The word "TECHNE" is written in a bold, sans-serif font and is enclosed within a horizontal oval border. This logo is centered on a thin horizontal line that spans the width of the page.

TECHNE

TC-3000 & TC-3000X

OPERATOR'S MANUAL

CONTENTS

	PAGE
INTRODUCTION	1
SAFETY AND INSTALLATION	2
SÉCURITÉ ET CONSIGNES D'INSTALLATION	5
SICHERHEITS- UND INSTALLATIONSINFORMATIONEN	8
INFORMAZIONI SULLA SICUREZZA E L'INSTALLAZIONE	11
INFORMACIÓN DE SEGURIDAD E INSTALACIÓN	14
BEFORE SWITCHING ON	17
CONTACT INFORMATION	18
TC-3000 & TC-3000X SPECIFICATIONS	19
USER OPERATION AND CONSUMABLES	21
Heated lid	21
Tubes or reaction vessels	22
0.5ml tubes	23
Switching on	23
FRONT PANEL CONTROLS	24
The LCD display	24
Indicators	24
The keys	25
PROGRAMMING THE TC-3000 & TC-3000X	26
Programming Options	27
Option 1: Creating programs using templates	27
Option 2: Creating a new program	31
Example programs	32
Programming	32
To program increment or decrement temperature or time.	35
To run a program	39
INFORMATION	41
CONFIGURATION	41
RUNNING THE TC-3000 & TC-3000X FROM A PC	44
Connecting one unit to the computer using an RS 232 cable	44
Connecting several units to the computer using the daisy chain cables	45
NOTES ON TRANSFERRING PROTOCOLS FROM OTHER THERMAL CYCLERS	47
USER MAINTENANCE	48
Cleaning your TC-3000/X	48
Heater Block removal	48
SOME FREQUENTLY ASKED QUESTIONS	49
ADDITIONAL INFORMATION	50
Fault Finding	50
Fuses	50
The heated lid over-temperature cut-out	50
Insulation testing	50
Interchangeable blocks	50
Accessories	51
Replacement parts	52

INTRODUCTION

The TC-3000 and larger capacity TC-3000X provide the researcher with the means of accurately controlling the temperature profile of samples. It has many scientific applications, including PCR¹ and sequencing. Both units can cycle samples between 4°C and 99°C.

The TC-3000 and TC-3000X are programmed by means of an integral keypad and LCD display. A program, which can be recalled from memory, consists of:

- a series of specified temperatures in °C
- the times for which each specified temperature will be held (Hold Times)
- the desired heating or cooling rates, in °C/sec, between each specified temperature
- optional instructions as to whether the times and/or the temperatures are to increase or decrease when a cycle is repeated.

The memory can store up to 80 programs.

Before using the thermal cycler, make sure you have read this manual carefully. If there is any doubt relating to the proper use of this equipment, the staff at Bibby Scientific Ltd or your supplier will be happy to assist you.

¹Purchase of this instrument conveys a limited non-transferable immunity from suit for the purchaser's own internal research and development and applied fields other than human in-vitro diagnostics under non-real-time thermal cycler patents of Applied Biosystems LLC.

SAFETY AND INSTALLATION

OPERATOR SAFETY

Please read this manual carefully before using the thermal cycler. If the equipment is not used in the manner described in this manual the protection provided by the equipment might be impaired.

All Bibby Scientific Ltd instruments are designed to conform to international safety requirements and are fitted with an over-temperature cut-out. If a safety problem should be encountered then switch off the unit at the mains socket and remove the plug from the electricity supply.

WARNING

HIGH TEMPERATURES ARE DANGEROUS as they can cause serious burns to operators and ignite combustible material. Users should be aware of the following potential hazards:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS
- DO NOT use combustible substances near hot objects
- DO NOT operate the instrument in the vicinity of inflammable liquids or gases
- DO NOT place any liquid directly into the instrument, always use the appropriate plastic consumables

After heating samples, remember that parts of the instrument, namely the lid and block and any associated accessories may be very hot.

UNPACKING

When unpacking please ensure that the following have been removed from the packaging:

- Thermal cycler
- Mains cable
- Guarantee card
- Decontamination certificate

The user is advised to keep the original packaging in case the instrument ever needs to be returned for service or repair. Bibby Scientific Ltd accepts no responsibility for damage incurred unless the unit is correctly packed and transported in its original packaging.

Please note that the screen of the thermal cycler is easily damaged by sharp objects such as pens, pencils and fingernails. This type of damage will be considered as misuse and invalidate the guarantee for this component.

INSTALLATION

The instrument should be carried using both hands. Never move or carry the instrument when in use or connected to the mains electricity supply.

1. All Bibby Scientific Ltd instruments are supplied with a power cable; this may be integral or plug-in.
2. Before connecting the instrument to the mains electricity supply, check the voltage against the rating plate (located on the back of the unit). Ensure that the voltage selector switch (located above

the On/Off switch) is set to the appropriate voltage for the local supply. Please note that the unit must be earthed to ensure proper electrical safety. (The TC-3000X has a switch-mode power supply which is suitable for use with universal AC input voltages between 90 and 260V. It is not necessary to change the fuses when using the TC-3000X). Connect the mains cable to a suitable plug according to the table below.

Connections	220V-240V, 50/60Hz Supply	110V-120V, 50/60Hz Supply
Live	brown	black
Neutral	blue	white
Earth	green/yellow	green

- UK ONLY: The fused plug supplied with the mains cable is fitted with a 10 Amp fuse to protect the instrument and the user.
- TC3000 units showing 230V, 50/60Hz on the rating plate also operate between 210 and 260V; units with 115V, 50/60Hz also operate between 110 and 130V and units marked with 100V will operate between 90 and 110V. In all cases the heating rate will degrade by approximately 8% at the extremes of the voltage range.
- Plug the mains cable into the socket on the back of the instrument.
- Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction. Position the instrument with a minimum distance all round of 200mm from walls or other items and between each unit if multiple units are being used to permit efficient air flow for each instrument.
- Switch on the instrument:
 - I mains switch On
 - O mains switch Off

Replacement cable

Should the mains lead need replacement, a cable of 1mm² of harmonized code H05VV-F connected to an IEC320 plug should be used. **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN.**

WORKING CONDITIONS

The thermal cyclers are designed to operate under the following conditions:

- Indoor use
- Ambient temperature range +5°C to +40°C
- Altitude to 2000m
- Mains supply fluctuations not exceeding 10%
- Over voltage category II IEC 60364-4-443
- Pollution degree 2

Note: The control specifications are quoted at an ambient temperature of 20°C. The specification may deteriorate outside an ambient temperature of 10°C to 30°C.

The instrument has been tested for radio frequency interference and is certified under EN61326.

GUARANTEE

The instrument is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is effective from the date of purchase; within this period all defective parts will be replaced free of charge provided that the defect is not the result of an accident, misuse or negligence. The screen of the thermal cycler can be damaged by sharp objects such as pens, pencils and fingernails. Damage of this sort will be considered misuse and will invalidate the guarantee for this component.

Servicing under this guarantee should be obtained from the supplier of the instrument.

This manual has been prepared for the convenience of Bibby Scientific Ltd's customers and nothing in this manual shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the unit or components.

Notwithstanding the description and specification(s) of the instruments contained in the operator's manual, Bibby Scientific Ltd reserves the right to make such changes as it sees fit to the instruments or to any of the components.

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
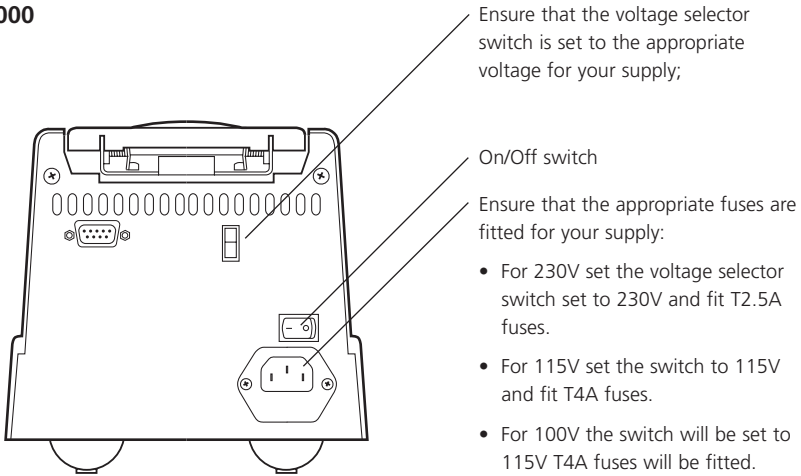
TC-3000X

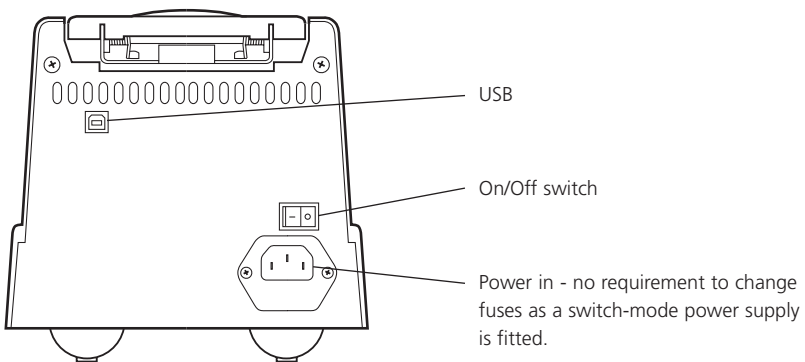CONTACT INFORMATION

For technical, sales or servicing information, contact your local Bibby Scientific Ltd dealer or:

UK and Rest of World

Bibby Scientific Ltd
Beacon Road, Stone
Staffordshire ST15 0SA, UK
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 813748
e-mail: techne@bibby-scientific.com
www.bibby-scientific.com

France

Bibby Scientific France SAS
ZI du Rocher Vert - BP 79
77793 Nemours Cedex
France
Tel: +33 1 64 45 13 13
Fax: +33 1 64 45 13 00
e-mail: bsf@bibby-scientific.fr
www.bibby-scientific.com

Middle East

Techne Middle East Ltd.
PO Box 27887, Engomi 2433
Nicosia
Cyprus
Tel: +357 22 660427
Fax: +357 22 660356
e-mail: sales@techneme.com

North and South America

Techne Inc.
3 Terri Lane, Suite 10
Burlington, N.J. 08016, USA
Tel: 800-225-9243
Fax: 609-589-2571
e-mail: labproducts@techneusa.com
www.techneusa.com

Italy

Bibby Scientific Italia Srl
Via Alcide de Gasperi 56
20077 Riozzo di Cerro al Lambro
Milano, Italia
Tel: +39 (0)2 98230679
Fax: +39 (0)2 98230211
e-mail: marketing@bibby-scientific.it
www.bibby-scientific.it

Far East

Bibby Scientific (Asia) Ltd
Room 607 Yen Sheng Centre
64 Hoi Yuen Road Kwun Tong
Kowloon Hong Kong
Tel: +852 3583 1581
Fax: +852 3583 1580
e-mail: bibby@bibby-scientificasia.com
www.bibby-scientificasia.com

TC-3000, TC-3000X SPECIFICATIONS

	TC-3000	TC-3000X
Temperature		
Temperature range	4°C to 99°C	4°C to 99°C
Temperature set point precision	0.1°C	0.1°C
Block uniformity (over full range)	± 0.1°C	± 0.2°C
Temperature accuracy	± 0.2°C	± 0.2°C
End of program cooling below ambient	Yes	Yes
Heating/Cooling Rate		
Heating rate, between 55°C and 90°C on a 50°C to 95°C step:		
0.2 ml micro tubes	3.6°C/sec	3.3°C/sec
0.5 ml micro tubes	3.6°C/sec	3.3°C/sec
Cooling rate, between 90°C and 55°C on a 95°C to 50°C step:		
0.2 ml micro tubes	2.0°C/sec	2.0°C/sec
0.5 ml micro tubes	1.8°C/sec	2.0°C/sec
Heated Lid		
Adjustable heated lid	No	Yes
Heated lid enable/disable	Yes	Yes
Selectable heated lid temperature	100°C to 115°C	100°C to 115°C
Over-temperature cut-out	Fixed at 145°C	Fixed at 145°C
Temperature sensor	Thermistor	Thermistor
Heater type	Etched foil	Etched foil
Heater power	33W	33W
Warm up time	4 min approximately (ambient to 105°C)	4 min approximately (ambient to 105°C)
The heated lid will only come on if the set temperature is above 35°C		
Programming		
Program naming using alpha keys	Yes	Yes
Program password protection	Yes	Yes
Number of programs (3 step)	80	80
Maximum number of steps per stage	35	35
Maximum number of stages per program	25	25
Minimum hold time	1 sec	1 sec
Maximum hold time	18 hrs	18 hrs
Incremental/decremental temperature	Yes	Yes
Incremental/decremental step hold time	Yes	Yes
Programmable ramp rate (heat or cool)	Variable in 0.1°C/sec steps	Variable in 0.1°C/sec steps
Programmable ramp rate resolution	0.1°C	0.1°C

	TC-3000	TC-3000X
Running Programs		
Pause button	Yes	Yes
Stop button	Yes	Yes
End of program alarm	Yes	Yes
Auto restart on power fail	Yes	Yes

Interconnectivity		
Connection to PC control program	Yes	Yes
Software updates from Techne web-site	Yes	Yes

Dimensions		
Height	185mm	185mm
Width	185mm	185mm
Length	335mm	335mm

Communication Ports		
PC connection to run "Gensoft" the PC control software.	RS232 serial	USB

Power		
Power consumption	230W	230W

The TC-3000 units are voltage selectable by a slide switch. When the unit is set to 115V it will work at any voltage between 110V and 130V. When the unit is set to 230V it will work at any voltage between 210V and 260V. Units marked 100V will work at any voltage between 90 and 110V.

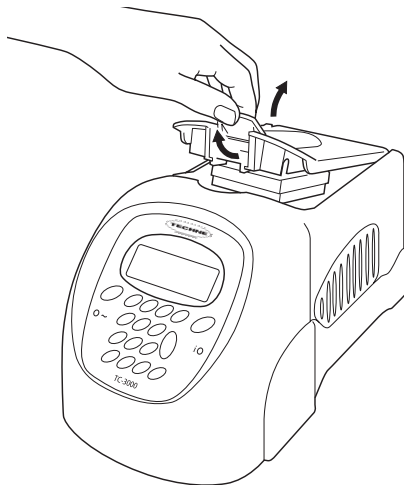
The performance will vary and will not necessarily meet the above typical specification at the extremes of voltage.

The TC-3000X unit is supplied with a switched mode power supplier so the unit automatically recognises the input voltage.

USER OPERATION AND CONSUMABLES

HEATED LID: TC-3000

To release the heated lid, lift the lid catch and the lid will open. To close the lid, push down gently until the catch clicks in place. **Do not close the lid without tubes being fitted in the block.**

HEATED LID: TC-3000X

The heated lid has a rotating knob to adjust the lid pressure on the samples, allowing for a variety of consumables to be used. To adjust the pressure:

- Adjust the orange rotating knob anticlockwise to give the least pressure.
- Place the samples in the block and place the lid in the down position.
- Gently rotate the knob clockwise until you can feel pressure beginning to be applied.
- Rotate the knob a further quarter of a turn and this will give the correct pressure.

Do not over-tighten the knob as this will cause deformation of the silicone rubber mat (where used) and reduce the effectiveness of the lid. Once the pressure has been set it does not need to be adjusted unless a different block or consumable is used.

The silicone rubber mat supplied with the unit (Part number 6106196) is for use with 48 well microplates when sealing using a plastic adhesive film. The mat should be placed on top of the sealing film to ensure good thermal contact with the heated lid.

TUBES OR REACTION VESSELS

Bibby Scientific Ltd does not recommend any specific tube or reaction vessel other than those described in this manual. We recommend using reaction volumes between 20 and 200 μ l. The tubes must withstand a pressure of 1 atmosphere at 100°C. Any vessel must be able to withstand the temperatures you are using without any danger of them deforming to the point where they fracture.

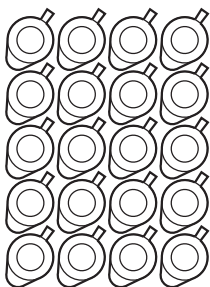
To test your tubes, put 25 μ l of water into each of 5 tubes and subject them to a typical thermal cycling protocol. At the end of the program, measure the volume remaining using a micropipette. A loss of more than 1 or 2 μ l indicates a vapour leak.

The amount of volume loss you observe and the change in reactant concentrations you can tolerate determine the minimum volume that can be used. Typical volume losses of 1 μ l in 30 cycles allow the use of samples of 20 μ l or less.

During the final cool-down, a ring of condensation may form above the liquid level but below the top of the sample block. This is normal and does not effect the reaction.

0.5ML TUBES

The heated lid mechanism of the TC-3000 with the 0.5ml block is optimised for flat top tubes; it is not suitable for use with domed cap tubes. In order to fit the maximum number of tubes in the block, they need to be placed so that all the hinges are in the same relative orientation. An example is illustrated below.

Place the tubes in the block in the same orientation in order to achieve the maximum capacity.

Always make sure the pressure on the heated lid is balanced by spreading the tubes across the block or inserting some dummy tubes if necessary.

PLATES AND SEALS

The 48 x 0.2ml block of the TC-3000X will accept 48-well plates as well as individual and strip 0.2ml tubes. To avoid the possibility of sample evaporation Bibby Scientific Ltd recommends that plates are sealed with appropriate caps or heat seals. We do not recommend the use of adhesive seals, but if their use is unavoidable then the supplied silicone rubber mat should be placed on top of the sealing film to ensure good thermal contact with the heated lid.

SWITCHING ON

TC-3000	
Version	x.xx
Block	25x0.2ml

When you switch on the unit a screen similar to the one shown here will appear briefly. This indicates the software version and block type installed in the unit.


Programs	[0]
Information	
Configuration	
[↑ ↓ , ← to select]	

When the initial screen has disappeared the screen shown here will automatically appear. There are three options:


1. 'Programs' for creating and running programs;
2. 'Information' for obtaining information about the unit e.g. serial number;
3. 'Configuration' for setting up program defaults and settings.

Each of these is explained in detail in the following sections.

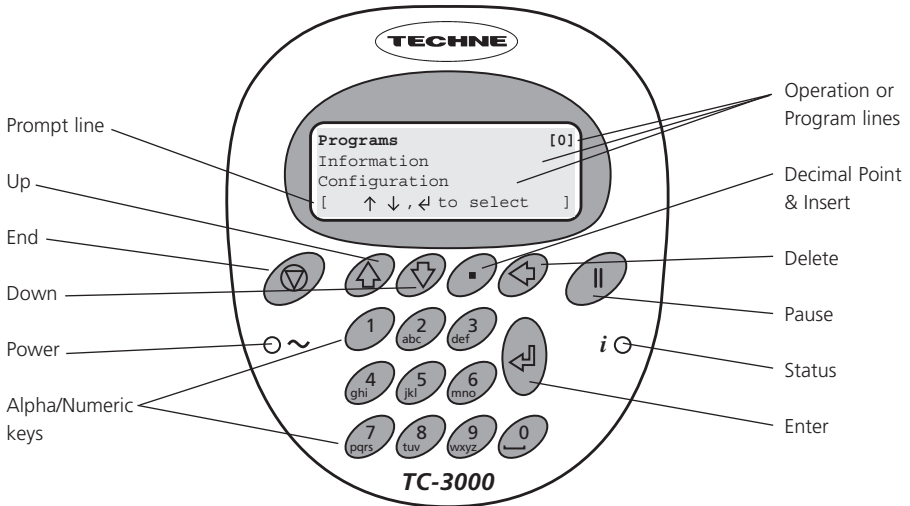
FRONT PANEL CONTROLS

THE LCD DISPLAY

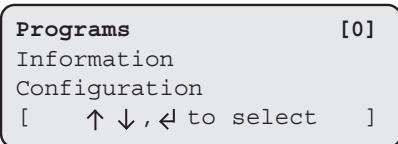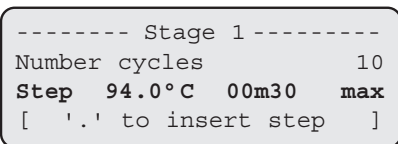The TC-3000 & TC-3000X have a four line LCD display. The top three lines are information or program lines, while the bottom line is a prompt line. The active line flashes on screen and is shown in bold in this manual.

[↑ ↓, ← to edit]

[← to delete step]

The bottom line may change or flash between two or three different prompts depending upon which program line is active. An example of a prompt line which flashes between three different prompts is shown here.

INDICATORS

○ ~ Power: This indicator is lit when there is power to the TC-3000 or TC-3000X.

i ○ Status: This indicator is lit when the unit is running a program. It is off when the unit is stopped. It will flash and beep slowly when the unit is paused. It will flash after a five second quick sequence of beeps when the unit reaches its final hold temperature.

THE KEYS

The End, Finish or Exit key

The "End" or "Finish" key is used to finish editing programs or to exit a program. It is also used to exit a sequence of key operations.

The Up arrow key

The "Up arrow" key is used for scrolling up through the lines on the screen. It is also used to increase certain fields when editing.

The Down arrow key

The "Down arrow" key is used for scrolling down through the lines on the screen. It is also used to decrease certain fields when editing.

The Decimal Point and Insert key

The "Decimal Point" key is used when editing numeric fields. It is also used as an Insert key during program editing to insert additional steps and other field sets.

The Delete key

The "Delete" key is used during program editing to delete steps and stages.

The Pause key

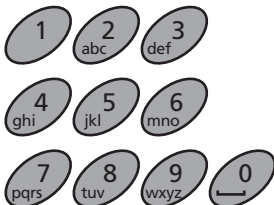
The "Pause" key is used during program editing to insert a pause after a stage or to insert increment and decrement times and temperatures. It is also used to pause a program when it is running.

The Enter or Accept key

The "Enter" key is used to enter an operation or the edit mode. It is also used to 'Accept' a field or program change.

The Alpha/Numeric keys

The "Alpha/Numeric" keys are used to enter numbers and letters during programming. If it is a numeric field then numbers will be inserted; if it is an alpha field then letters will be inserted. Each key will cycle through the letters shown on the key (0 will insert a space). To insert a number, press and hold the appropriate key.

PROGRAMMING THE TC-3000 & TC-3000X

The table below illustrates the symbols which appear on the LCD display and the relevant keypad button to which they refer.

	In the prompt line is equivalent to the "End" key.	
	On a program line means the program is locked; it can be copied and then edited but not edited directly.	
	Up arrow	
	Down arrow	
	Enter	
	Pause	

The following terms are used when programming the thermal cyclers:

- Stage** A 'Stage' is the name given to the main repeated part of a program containing the 'Steps' to be repeated. A program can contain a maximum of 25 stages although typically only one or two are used.
- Cycle** A 'Cycle' is the number of repetitions of a 'Stage'
- Step** A 'Step' is a user defined program line containing a temperature, a hold time and ramp rate. Typically three steps are created within one 'Stage'.

```

Programs                                [0]
Information
Configuration
[   ↑ ↓ , ← to select   ]
    
```

To enter the **programming** section of the software, highlight 'Programs' using the "Up Arrow" or "Down Arrow" keys ↑ ↓

Press "Enter" ←

```

* * New Program * * [ ]
2 STEP TEMPLATE  501  9
3 STEP TEMPLATE  502  9
[   ↑ ↓ , ← to select   ]
    
```

This screen lists all the programs saved on the unit. Use the "Up Arrow" and "Down Arrow" keys to scroll through the list. ↑ ↓

The number next to the program name is the unique "Speed dial" number allocated to the program for rapid retrieval (see below).

PROGRAMMING OPTIONS

There are two options available for creating programs on the TC-3000 & TC-3000X:

1. Copy and edit one of the pre-written templates.
2. Create a new program from the defaults.

OPTION 1: CREATING PROGRAMS USING TEMPLATES

There are two pre-written basic PCR templates: a two-step and a three-step template. These can be copied and modified as required.

2 STEP TEMPLATE

```

Heated Lid          105°C
Preheat Lid         on
Pause              off
-----
In denat            94°C 5m00
Hot Start           off
----- Stage 1-----
Number cycles       30
Step 94.0°C 0m30 max
Step 68.0°C 1m00 max
-----
Final Extn 72.0°C 5m00
Final hold          10.0°C
    
```

3 STEP TEMPLATE

```

Heated Lid          105°C
Preheat Lid         on
Pause              off
-----
In denat            94.°C 5m00
Hot Start           off
----- Stage 1-----
Number cycles       30
Step 94.0°C 0m30 max
Step 55.0°C 0m30 max
Step 72.0°C 0m30 max
-----
Final Extn 72.0°C 5m00
Final hold          10.0°C
    
```

Other templates available include Ice bucket (incubation at 4°C), Ligation (incubation at 15°C) and an RT-PCR template.

```

* * * New Program * * *
2 STEP TEMPLATE 501
3 STEP TEMPLATE 502
[ ↑ ↓, ← to select ]
    
```

Use the "Down Arrow" key ↓ to select an existing 'Template' program from the menu, e.g. '2 Step Template' or '3 Step Template'.

Press the "Enter" key ↵ to accept it.

```

2 STEP TEMPLATE [501]
Run program
View program
[ ↑ ↓, ← to select ]
    
```

The Template is already defined and can be run as it is or easily copied, modified and saved under a new name.

```

Run program
View program
Copy program
[ ↑ ↓, ← to select ]
    
```

Use the "Down Arrow" key ↓ to select 'Copy Program' Press the "Enter" key ↵ to accept it.

```

      UNNAMED #
Heated Lid      105°C
Preheat Lid      on
[   ↑ ↓ , ⏮ to edit   ]
    
```

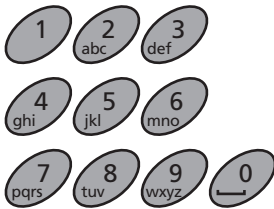
A new name can now be given to the copied program (see below).

To give a program a name

```

■      UNNAMED #
Heated Lid      105°C
Preheat Lid      on
[   ↑ ↓ , ⏮ to edit   ]
    
```

To type in a name for the program, press "Enter" ⏮ and a cursor will appear in the position shown.

Using the keypad, type in the name you wish to give the program.

The letters will scroll round with successive presses; for example the "2" key will give 'A' with one press; 'B' with two presses; 'C' with three presses and '2' with four presses. Five presses will again give 'A'.

When the name is complete, press "Enter" ⏮

```

      DONNA
Heated Lid      105°C
Preheat Lid      on
[   ↑ ↓ , ⏮ to edit   ]

[   ▽ to save   ]
    
```

There are now two options:

1. Save the program and run it as it is or
2. Edit the program then save and run it.

To save the program

```

      DONNA
Heated Lid      105°C
Preheat Lid      on
[   SAVE?   ]

[   ⏮ YES or ▽ NO   ]
    
```

A program can only be run once it has been saved. Press the "End" ▽ key.

The prompt line will flash between 'SAVE?' and 'YES or NO'.

Press "Enter" ⏮ for 'Yes' to save the changes you have made.

If you press "End" ▽ for 'No' then all the changes you have made will be lost.

```

      DONNA      [7]
Run Program
View Program
[   ↑ ↓ , ⏮ to select   ]
    
```

The program is now ready to run and will be allocated with a unique 'Speed dial' number which can be used to rapidly locate the program from the main start-up screen.

"Up Arrow" or "Down Arrow" keys ↑ ↓ to scroll through the remaining list of options.

To edit a field or step in an existing program

```

Programs [0]
Information
Configuration
[  ↑ ↓ , ↵ to select  ]
    
```

```

DAVID 8 [D]
DONNA 7
[  ↑ ↓ , ↵ to select  ]
    
```

When the unit is switched on the display will show this screen.

If known, type in the 'Speed dial' number for the required program and press "Enter" ↵. This will take you directly to the required program.

If you do not know the 'Speed dial' number, press "Enter" ↵ to access the list of stored programs.

Either:

1. Scroll to the required program using the "Up Arrow" or "Down Arrow" keys ↑ ↓ or
2. Type the first letter of the program name in the 'New program' field; this will then list all the program names beginning with that letter. Use the "Up Arrow" or "Down Arrow" keys ↑ ↓ to select the required program.

Press "Enter" to accept ↵

```

DONNA [7]
Run program
View program
Edit program
Copy program
Delete program
Set password
[  ↑ ↓ , ↵ to select  ]
    
```

The full set of options for a program is shown here; these are shown on the LCD screen three lines at a time.

Use the "Down Arrow" and/or the "Up Arrow" keys ↑ ↓ to select the option you require.

Press the "Enter" key ↵ to select the option.

Edit a selected field or step

```

          DONNA
Heated Lid 105°C
Preheat Lid on
[  ↑ ↓ , ↵ to edit  ]
    
```

Select 'Edit program' and a screen similar to this will appear.

Use the "Down arrow" and/or the "Up arrow" keys ↑ ↓ to select the field or step you want to edit.

You can now edit the active field or step by first pressing the "Enter" key ↵

```

[  ▽ to save  ]
    
```

```

DONNA
Heated Lid      105°C
Preheat Lid      on
[   edit temperature   ]

[   ← to disable   ]

```

To edit a temperature, type in the required temperature. In the example here, the heated lid temperature is being edited.

Press "Enter" ↵ to accept.

Use the "Down Arrow" and/or the "Up Arrow" keys ↑ ↓ to select the next field or step you wish to edit.

```

DONNA
Heated Lid      105°C
Preheat Lid      on
[   ↑ ↓ to change   ]

[   turn on/off   ]

```

To edit a field with on/off settings, first select the field and press "Enter" ↵

Use the "Up Arrow" and "Down Arrow" keys to toggle ↑ ↓ between the settings.

Press "Enter" ↵ to accept.

```

DONNA
Preheat Lid      on
Sample cooling     off

```

Sample cooling is a new function on the TC-3000X that enables to be actively cooled to 4°C whilst the heated lid is reaching the correct temperature.

```

Pause      off
-----
In denat   94.0°C   00m30
[ edit temperature ]

[   ← to disable   ]

```

To edit a step with both a temperature and a time, first press "Enter" ↵ and type in the new temperature. Press "Enter" ↵ again. Now type in the new time. The default is minutes.

1. For whole minutes, type the number required then "Enter" ↵.
2. To toggle between minutes and hours, use the "Decimal Point" key.
3. For values less than 1 minute, first press the "Decimal Point" key followed by the number of seconds required.
4. For minutes and seconds, type the number of minutes followed by the "Decimal Point" key then the number of seconds.

```

Pause      off
-----
In denat   94.0°C   00m30
[ edit time   ]

[   .for hrs/mins   ]

```

Once the values have been added, press "Enter" ↵ to accept.

```

-----
Fin extn 72.0°C   5m00
Fin hold  10.0°C
[   ↑ ↓, ↵ to edit   ]

[   ▽ to save   ]

```

Once you have finished editing and you are happy with the changes, press the "End" key ▽ followed by the "Enter" key ↵ to save.

If you press the "End" key ▽ a second time for 'No' then all the changes you have made will be lost.

To copy a selected program

UNNAMED #	
Heated Lid	105°C
Preheat Lid	on
[↑ ↓ , ← to edit]	
[▽ to save]	

If you select 'Copy program' a screen similar to this will appear and the software will automatically give the program the next number until it is named.

Press "Enter"  and type in a new name. You will then be able to edit the program as previously described.

To delete a selected program

DONNA	
delete program	
SURE?	
[← Yes or ▽ No]	

If you select 'Delete program' a screen similar to this will appear.

Press "Enter"  to delete the program or "End"  to return to the program screen.

To set a password for a selected program

DONNA	
New passwd	■
[← to accept]	
[▽ to exit]	

If you want to password protect the program so that other users cannot edit it, select 'Set password'. A screen similar to this will appear. Type in the password you wish to give it and press "Enter" . The program list will then display the key symbol  next to the name to show that it is locked. Other users will still be able to copy/view your program but not edit it.

OPTION 2: CREATING A NEW PROGRAM

A second option is to create a completely new program without using templates.

UNNAMED #	
Heated Lid	105°C
Preheat Lid	on
Pause	off

In denat 94.°C	5m00
Hot Start	off

Final Extn 72.0°C	5m00
Final hold	10.0°C

A number of typical values are pre-entered including:

Heated lid temperature
Preheat lid
Sample cooling (TC-3000 shown in diagram)
Initial denaturation
Hot start
Final extension
Final hold

These values can be edited and the additional stages and steps required for cycling can be added by following the instructions detailed on the following pages.

EXAMPLE PROGRAMS

A couple of example thermal cycling programs are shown below.

Typical 3-Step PCR

```

Heated Lid          105°C
Preheat Lid         on
Pause               off
-----
In denat 94.0°C    5min
Hot Start           off
----- Stage 1 -----
Number cycles       30
Step 94.0°C        0m30    max
Step 55.0°C        0m30    max
Step 72.0°C        0m30    max
-----
Final Extn         72.0°C    5m
Final hold         4°C      10m
  
```

3-Step PCR with multiple 'Stages'

```

Heated Lid          105°C
Preheat Lid         on
Pause               off
-----
In denat 94.0°C    5min
Hot Start           off
----- Stage 1 -----
Number cycles       10
Step 94.0°C        0m30    max
Step 60.0°C        0m30    max
Step 72.0°C        0m30    max
-----
Number cycles       20
Step 94.0°C        0m30    max
Step 55.0°C        0m30    max
Step 72.0°C        0m30    max
-----
Final Extn         72.0°C    5m
Final hold         4°C      10m
  
```

PROGRAMMING

```

* * New Program * * [ ]
2 STEP TEMPLATE 501  ⏴
3 STEP TEMPLATE 502  ⏴
[  ↑ ↓, ⏴ to select ]
  
```

```

          UNNAMED #
Heated Lid          105°C
Preheat Lid         on
[  ↑ ↓, ⏴ to edit ]

[  ▾ to save ]
  
```

Create a new program by selecting ** New Program ** from the menu and press "Enter" ⏴

Give the program a name and edit any of the default fields as described above.

To add a stage and then edit it

```

In denat 94.0°C    5m00
Hot Start           off
-----
[  ⏴ to insert stage ]
  
```

Use the "Down Arrow" ↓ key to scroll to the dotted line below 'Hot Start'; it will start to flash.

Press "Enter" ⏴ to insert a stage.

```

----- Stage 1 -----
Number cycles          █
Step  94.0°C 00m30    max
[    ␣ edit no.cycles  ]

```

A new stage will be inserted which consists of a 'Number cycles' field and one step.

Edit the number of cycles by typing the number required then press "Enter" ␣

```

----- Stage 1 -----
Number cycles          10
Step 94.0°C 00m30    max
[    '.' to insert step  ]

[    ↑ ↓, ␣ to edit      ]
[    ← to delete step    ]

```

Use the "Down Arrow" ↓ key to move to the step.

Press "Enter" ␣ to edit the step as described above.

Press "Enter" ␣ again to accept.

To insert additional steps

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C 00m30    max
[    '.' to insert step  ]

[    ↑ ↓, ␣ to edit      ]

```

To insert a new step before the first step, use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to select 'Number cycles'.

Press the "Decimal Point" key to insert a new step.

To insert a new step after an existing step, use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to select the step.

Press the "Decimal Point" key to insert a new step.

To delete a step

```

----- Stage 1 -----
Number cycles          10
Step 94.0°C 00m30    max
[    '.' to insert step  ]

[    ↑ ↓, ␣ to edit      ]
[    ← to delete step    ]

```

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to select the step you wish to delete.

Press "Delete" ← to delete the step.

```

----- Stage 1 -----
Number cycles          10
Step 94.0°C 00m30    max
[    DEL STEP ?         ]

[    ␣ Yes or ▽ No      ]

```

The prompt line will flash between 'DEL STEP?' and 'YES or NO'.

Press "Enter" ␣ to delete the step.

or

Press "End" ▽ to return to editing the step.

To add a new stage before an existing one

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C  00m30  max
[    ⏮ to insert stage ]

[    ⏪ to delete stage ]

```

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to highlight the stage number.

Press "Enter" ⏵

A new stage will be added before the existing stage. The other stage numbers will be adjusted as appropriate. The new stage can now be edited as described above.

To add a new stage at the end

```

-----
Final Extn 72.0°C      5m00
Final hold           10.0°C
[    ⏮ to insert stage ]

```

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to highlight the dotted line below the last stage.

Press "Enter" ⏵

A new stage will be added with the next consecutive number. The new stage can now be edited as described above.

To delete a stage

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C  00m30  max
[    ⏮ to insert stage ]

[    ⏪ to delete stage ]

```

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to highlight the stage you wish to delete.

Press "Delete" ⏴

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C  00m30  max
[    DEL STAGE?      ]

[    ⏮ Yes or ▽ No   ]

```

The prompt line will flash between 'DEL STAGE?' and 'YES' or 'NO'.

Press "Enter" ⏵ to delete the stage.

To insert a 'pause' into a program

```

Step  94.0°C  00m30  max
----- Stage 2 -----
Number cycles          █
Step  94.0°C  00m30  max

[  '.' to insert step  ]
[  ↑ ↓ , ← to edit   ]
    
```

To insert a pause after an existing stage, follow the steps for inserting a stage as above.

Highlight the 'Number cycles' line then press "Pause" | |

```

Step  94.0°C  00m30  max
----- Stage 2 -----
Number cycles          0
Step  94.0°C  00m30  max

[  CHANGE TO PAUSE    ]
[  ← Yes or ▽ No     ]
    
```

The prompt line will flash between 'CHANGE TO PAUSE?' and 'YES or NO'.

Press "Enter" ← to insert a pause.

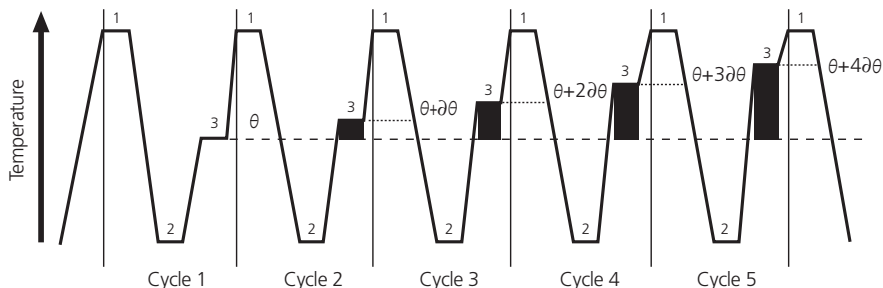
```

----- Stage 2 -----
Pause
-----
[  ← to insert step  ]

[  ← to delete step  ]
    
```

The stage number will then be highlighted with 'Pause' as the only line.

TO PROGRAM INCREMENT OR DECREMENT TEMPERATURE OR TIME


Under normal circumstances, the Hold Temperature of all steps is constant. However, it is possible to automatically increment or decrement the temperature of a specified step of a programme.

Temperature profile of incremented Hold Temperature at step 3 of each cycle

The Hold Temperature of the incremented/decremented step is the defined Hold Temperature plus the summation of the increments/decrements. The first cycle is never incremented/decremented, only subsequent cycles. If you select a decrement, the Hold Temperature is prevented from falling below 4°C. If you select an increment, the Hold Temperature is prevented from rising above 99°C.

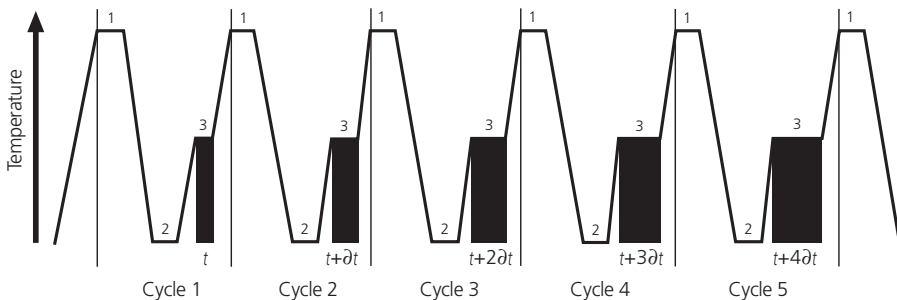Similarly, under normal circumstances, the Hold Time of all steps is constant. However, it is possible to automatically increment or decrement the duration of a specified step of a program.

Temperature profile of incremented Hold Time at step 3 of each cycle

The Hold Time of the incremented/decremented step is the defined Hold Time plus the summation of the increments/decrements. The first cycle is never incremented/decremented, only subsequent cycles. If you select a decrement, the Hold Time is prevented from falling below one second.

Number cycles	30
Step 94.0°C 00m30	max
Step 58.0°C 00m30	max
[, to edit]	

[← to delete step]
[`.' to insert step]

Number cycles	30
Step 94.0°C 00m30	max
Step 58.0°C 00m30	max
[CHANGE STEP ?]	

[← Yes or ▽ No]

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to move to the step where you want an increment or decrement temperature or time

Press "Enter" ↵ to go into edit mode.

Press "Pause" | |

The prompt line will change to flash between 'CHANGE STEP?' and 'YES' or 'NO'.

Press "Enter" ↵

The cursor will again move into the edit mode.

Now press the "Up Arrow" key ↑

```

Number cycles          30
Step  94.0°C   00m30   max
  fst  55.0°C   00m30   max
[    ↑ ↓, ← to edit  ]
  
```

[▽ to save]

```

Number cycles          30
Step  94.0°C   00m30   max
  fst  65.0°C   00m30   max
[    edit temperature  ]
  
```

```

Step  94.0°C   00m30   max
fst   65.0°C   00m30   max
lst   55.0°C   00m30
[    ↑ ↓, ← to edit  ]
  
```

[← to delete step]

[`.' to insert step]

```

Step  94.0°C   00m30   max
fst   65.0°C   00m30   max
lst   55.0°C   1m30
[    edit time        ]
  
```

[.For hrs/mins]

The step will change to 'fst', meaning "first", and the software will insert another line after it entitled 'lst', meaning "last".

Press "Enter" ↵ twice to go into edit mode.

Use the "Alpha/numeric" keys to change the values of the 'fst' line to the first temperatures and/or times you require.

Press "Enter" ↵ after each field you change.

You will exit 'Edit Mode' when you reach the end of the line.

Use the "Down Arrow" key ↓ to move onto the 'lst' line.

Press "Enter" ↵ to edit the step.

Use the "Alpha/numeric" keys to change the values of the 'lst' line to the final temperatures and/or times you require.

Press "Enter" ↵ for each field you change.

You will exit 'Edit Mode' when you reach the end of the line.

When you run the program it will automatically increment or decrement in equal steps from the first time and/or temperature to the last over the number of cycles in the stage. For example: if the first temperature is 65°C and the last temperature is 55°C, over 30 cycles the temperature will decrease 0.33°C each cycle. Similarly, in the example shown, the time would increase from 30 seconds to 1 minute 30 seconds in 2 second increments.

To save the program

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C  00m30  max
[      SAVE?      ]

[    ↵ Yes or ▽ No    ]
    
```

A program can only be run once it has been saved. Press the "End" ▽ key.

The prompt line will flash between 'SAVE?' and 'YES or NO'.

Press "Enter" ↵ for 'Yes' to save the changes you have made.

If you press "End" ▽ for 'No' then all the changes you have made will be lost.

The program is now ready to run and will be allocated with a unique "Speed dial" number which can be used to rapidly locate the program from the main start-up screen.

Use the "Up Arrow" and "Down Arrow" ↑ ↓ keys to scroll through the list of options.

```

                GEORGE      (10)
Run Program
View Program
[  ↑ ↓ , ↵ to select  ]
    
```

Stopping what you are doing

At any time during programming the TC-3000 or TC-3000X, when you are satisfied with the program, you can finish and accept what you have done by first pressing the "End" ▽ key.

Then follow the instructions on screen. If you are not sure, wait for the instructions in the prompt line to scroll back to the first line that you see.

TO RUN A PROGRAM

```

Programs [0]
Information
Configuration
[ ↑ ↓ , ↵ to select ]
    
```

```

Programs [7]
Information
Configuration
[ ↑ ↓ , ↵ to select ]
    
```

```

  * * New Program * * [ ]
2 STEP TEMPLATE 501  3
3 STEP TEMPLATE 502  3
[ ↑ ↓ , ↵ to select ]
    
```

```

DAVID 8 [D]
DONNA 7
[ ↑ ↓ , ↵ to select ]
    
```

```

DONNA [7]
Run Program
View Program
[ ↑ ↓ , ↵ to select ]
    
```

```

DONNA

[ RUN PROGRAM? ]

[ ↵ Yes or ▽ No ]
    
```

When the unit is switched on the display will show this screen.

If known, type in the 'Speed dial' number for the required program and press "Enter" ↵. This will take you directly to the required program.

If you do not know the 'Speed dial' number, press "Enter" ↵ to access the list of stored programs.

Next either:

1. Scroll down to the required program using the "Up Arrow" and "Down Arrow" ↑ ↓ keys.
2. Or type the first letter of the program name in the "New program" field; this will then list all the program names beginning with that letter. Use the "Up Arrow" and "Down Arrow" keys ↑ ↓ to select the required program.

Press "Enter" to accept ↵ For example 'Donna'.

To run the program, select 'Run program' and press "Enter" ↵

A screen similar to this one will appear.

Press "Enter" ↵ to start the run.

Cycle 4/30	56m08
Sample 58.0°C	
Hold @ 58.0°C	00m17
[▽ to STOP]	

[|| to PAUSE]
[↵ for info]

During the run the LCD screen will report:

1. The cycle number
2. The time remaining (counting down in seconds)
3. The sample temperature (whether holding or ramping)
4. The current step of the run
5. A count down of the hold time.

An example is shown here.

Manual pause

Cycle 24/30	19m26
Sample 72.0°C	
Hold @ 72.0°C	00m10
[PAUSE?]	

[↵ Yes or ▽ No]

To pause a program while it is running, press "Pause" || The run will continue but the prompt line will alternate 'PAUSE?' and 'YES or NO'

To confirm that you want the program to pause press "Enter" ↵

If you have pressed the "Pause" key by mistake and you want the program to continue, press "End" ▽

The unit will emit a beeping sound while paused.

To resume the run, press "Pause" || followed by "Enter" ↵ to continue.

Cycle 24/30	19m26
Sample 72.0°C	
Paused	
[▽ to STOP]	

[|| to CONTINUE]
[↵ for info]

Manual stop

Cycle 24/30	10m39
Sample 94.0°C	
Hold @ 94.0°C	00m10
[STOP PROGRAM?]	

[↵ Yes or ▽ No]

To stop a program while it is running, press "End" ▽

The run will continue but the prompt line will alternate 'STOP PROGRAM?' and 'YES or NO'

To confirm that you want to stop the program press "Enter" ↵

If you have pressed the "End" ▽ key by mistake and you want the program to continue, press "End" ▽ again.

When a program is stopped the unit will emit a beeping sound and the original program screen will appear.

Program Information

```

DONNA
Heated Lid           on
Status               ramping
[      ↵ to exit    ]

```

Pressing "Enter" ↵ while a program is running will give information about the current run including which program is actually running and the unit status e.g. holding or ramping.

An example is shown here.

INFORMATION

```

Programs              [0]
Information
Configuration
[  ↑ ↓, ↵ to select  ]

```

The **Information** field on the front screen contains all the unit-specific information such as serial number and software version.

Press "Enter" ↵ to access the unit information.

```

Block                20x0.5ml
Block s/n            139160-2
Unit s/n              GP0000-01
S/w version           14.zz
H/w version           01
Comms address         1
Unit cycles           624
Block cycles          69

```

The complete list is displayed here.

Note that the "Comms address" must be unique for each unit if more than one cycler is linked to a PC using the Gensoft control software.

CONFIGURATION

Warning! The Configuration software should only be accessed by designated personnel, as some parameters can be changed or may be lost when using it.

```

Programs              [0]
Information
Configuration
[  ↑ ↓, ↵ to select  ]

```

The **Configuration** field on the front screen allows a system administrator access to set up all the default settings for the programs, time and date settings and to change and remove passwords.

```

CONFIGURATION
Password             pad■
[      ↵ to accept    ]

[  ▽ to exit          ]

```

Access to the 'Configuration' set up is password protected.

To enter configuration, use the alpha keys to type in 'pad'

Press "Enter" ↵

```

Program defaults
Settings
Passwords
Controller reset

```

Program defaults

```

Program defaults
Settings
Passwords
Controller reset

```

```

Heated Lid           105°C
Preheat Lid           off
Pause                 off
-----
In denat      94°C    5m00
Hot Start                     off
-----
Final Extn 72.0°C    5m00
Final hold           10.0°C
[  ↑ ↓ , ← to select  ]

```

There are a number of options which can be changed; these are listed here.

Use the "Up Arrow" and "Down Arrow" ↑ ↓ keys to select then press "Enter" ↵ to access any of the options. Each option is discussed in detail below.

Highlight 'Program defaults' then press "Enter" ↵

The complete list of program defaults is given here. These are the values that all new programs will take when you start to create a new program.

To edit any line use the "Up Arrow" or "Down Arrow" keys ↑ ↓ to select the line you wish to change.

Press "Enter" ↵ then change the value using the numerical keys. Press "Enter" ↵ to accept the new value.

To edit a field with on/off settings, first select the field and press "Enter" ↵

Use the "Up Arrow" and "Down Arrow" ↑ ↓ keys to toggle between the settings. Press "Enter" ↵ to accept.

When you have set all the defaults, press the "Enter" ↵ key to go back to the Configuration menu.

Note: altering the default values does not change existing programs.

Settings

```

Program defaults
Settings
Passwords
Controller reset

```

Highlight 'Settings' then press "Enter" ↵

```

Set date      xx / xx / xx
Set time      xx : xx
Date format    dmy
Auto restart   on
Beeper        on
Coms address   1
Coms speed    4800
[  ↑ ↓ , ← to select  ]

```

The complete list of settings is given here.

Auto restart "on" will allow the unit to restart from where it left off should there be a power failure and the power returns.

The "Comms address" refers to the location of a unit in a daisy chain and must be unique if there are several units linked to each other.

The optimum "Coms speed" is set at 4800. When using the Gensoft control software, ensure all cyclers are set to the same value.

Edit any line as described for 'Program defaults' above. Press "Enter" ← to accept the new value.

When you have finished, press the "Enter" ← key to go back to the 'Configuration' menu.

Passwords

```

Program defaults
Settings
Passwords
Controller reset

```

The 'Passwords' function allows the administrator to change the 'Admin password' (which allows access to Configuration) and also to delete any passwords associated with programs.

Highlight 'Passwords' then press "Enter" ←

```

* * Admin passwd * *
DONNA
GEORGE
[  ↑ ↓ , ← to select  ]

```

The screen will appear as opposite.

We suggest that you change the 'Admin password' at this stage in order to protect the configuration settings.

```

      ADMIN PASSWD
Password  █
Confirm
[  ← to accept  ]

```

Use the "Up Arrow" ↑ key to highlight 'Admin passwd

'Press "Enter" ← and type in the new password. Repeat to confirm and accept.

The screen will go back to the 'Configuration' menu.

```

* * Admin passwd * *
DONNA
GEORGE
[  DEL PASSWORD?  ]
[  ← Yes or ▽ No  ]

```

To delete a password associated with a program, use the "Up Arrow" and "Down Arrow" ↑ ↓ keys to highlight the program.

Press "Enter" ←

Press "Enter" ← again to delete the password.

Controller reset

```

Program defaults
Settings
Passwords
Controller reset
    
```

The 'Controller reset' will delete all saved programs from the unit.

Highlight 'Controller reset' then press "Enter" ↵

```

          CONTROLLER RESET
This will delete
all programs
[          SURE ?          ]
[  ↵ Yes or ▽ No  ]
    
```

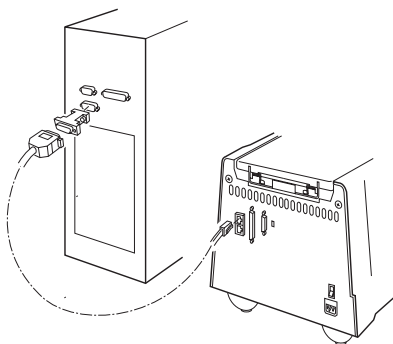
To delete all the programs, press "Enter" ↵

RUNNING THE TC-3000 & TC-3000X FROM A PC


The thermal cyclers can be run from a PC using Bibby Scientific Ltd's Gensoft software. This is available to download from the Techne web-site at www.techne.com. Instructions for using Gensoft are included in a separate manual.

Connecting one unit to the computer using an RS 232 cable

If you are driving one TC-3000 unit you will need part number FGEN232 which includes a 9-pin female to female, null modem, RS232, fully screened cable, Gensoft on disk and an instruction sheet. To connect a single TC-3000X unit, connect the unit to the PC using the USB cable, part number 6205046.

TC-3000

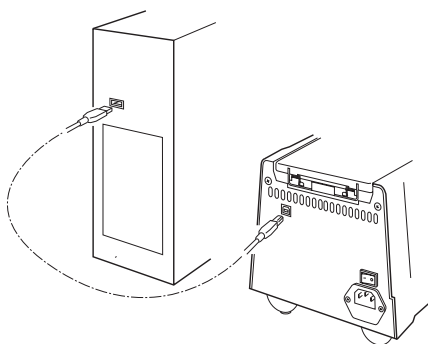
TC-3000X

Connecting several TC-3000 units to the computer using the daisy chain cables

If you are driving two or more units then you will need part number FGEN485D, a 230V UK power pack set, or FGEN485E, a 230V European power pack set, or FGEN485P, a 120/100V USA power pack set. A power pack set includes: the connection to a computer (which also converts to RS485), a power pack, the necessary cables, a termination box and the software on disk with an instruction sheet. You will also need one or more of the following:


-
- FGEN TWO** A two unit set which includes two connectors for the cyclers and a cable to interconnect them.
 - FGEN FIVE** A five unit set which includes five connectors for the cyclers and four cables to interconnect them.
 - FGEN TEN** A ten unit set which includes ten connectors for the cyclers and nine cables to interconnect them.
 - FGEN ONE** A one unit set for extending the daisy chain further which includes one connector for the cycler and one cable for connecting it to the daisy chain.
-

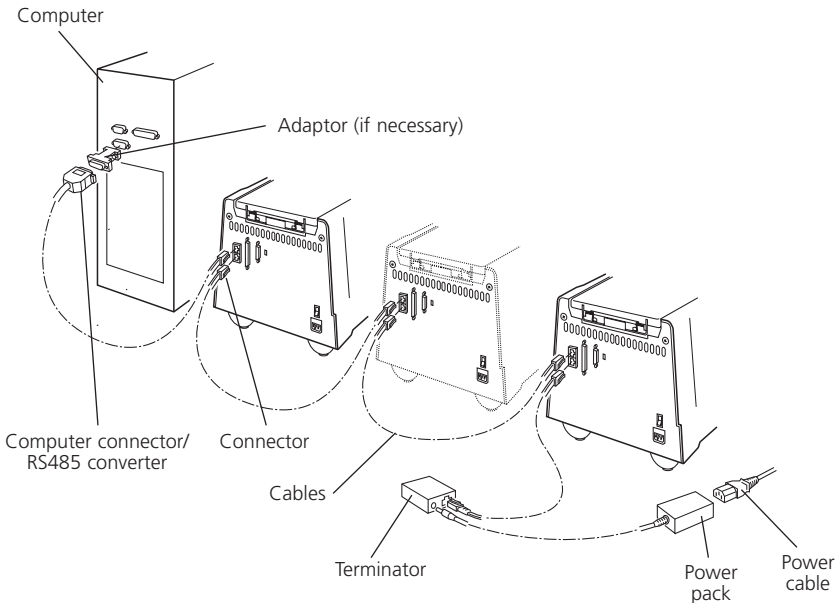Notice from the diagram that the cable connections to the cyclers are alternately bottom to top on the connector.

Connecting several TC-3000X units to the PC.

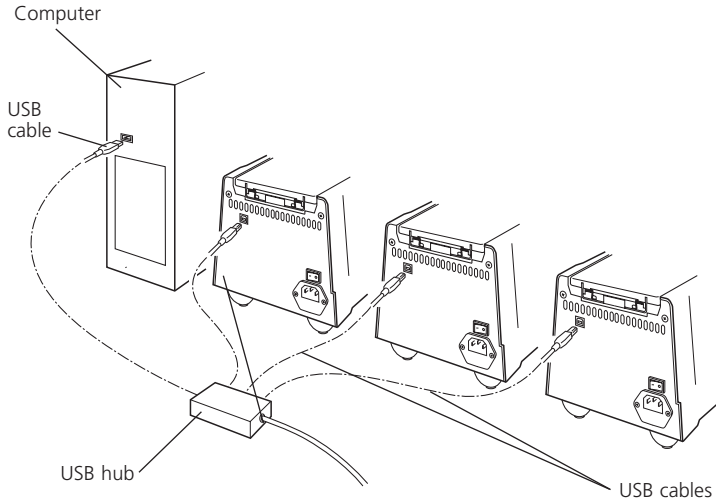If you are driving two or more units then you will need a USB cable for each unit, part number 6205046 plus a hub, which will connect multiple units together.

NOTES ON TRANSFERRING PROTOCOLS FROM OTHER THERMAL CYCLERS

The TC-3000 and TC-3000X can heat up to 3.6°C/s and provides the highest levels of temperature uniformity. Many slower machines have historically begun the countdown on hold times before the set temperature is reached in order to avoid any overshooting in sample temperature. The slower the heating rate the more significant the loss in hold time, i.e. if a 30 second hold time is set at 55°C then perhaps 5 to 10s will be lost in between the start of the time countdown and actually reaching the set temperature. The overall effect of this method of control is that your sample, if cooling from a higher temperature, is subjected to an environment with an elevated temperature compared to the target temperature for a significant part of the hold time.

The TC-3000 and TC-3000X provide the user with 'what-you-program-is what-you-get'. Due to the rapid and responsive nature of the Peltier heating system, the time between the timer countdown initiating and actually reaching the set temperature has been reduced to as little as 1s. This ensures that the sample is only subjected to the exact temperature required.

Taking the above into consideration it must therefore be noted that care must be taken when re-optimizing protocols from a slower machine.

Trouble Shooting

1. The optimum annealing temperature may in fact be higher than the set temperature used in the slower machine. Try increasing it by 1 or 2°C.
2. The hold time can be reduced once the optimum temperature has been found. Try reducing it by 5 to 10s.

USER MAINTENANCE

CLEANING YOUR TC-3000 OR TC-3000X

Before cleaning your unit, disconnect from the power supply and allow it to cool to below 50°C.

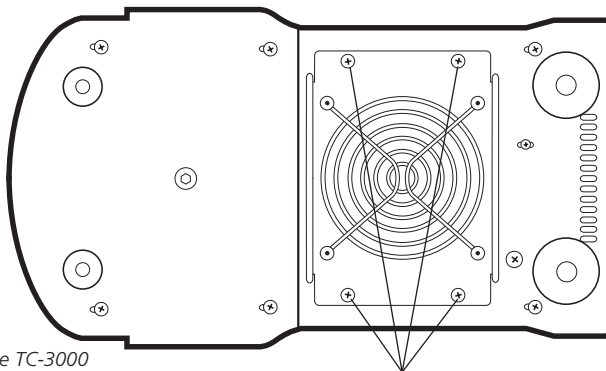
CLEANING YOUR BLOCK

The heating/cooling block, including wells and flat surfaces, should be cleaned regularly to ensure optimum heat transfer to the samples. Always clean the block if there has been a spillage. Use a cloth or cotton buds dipped in a fresh, 50:50 water/isopropanol solution and make sure that no deposits are left in the wells.

In the case of radioactive spillages the heater block can be removed from the unit for more detailed cleaning. Bibby Scientific Ltd recommends that you use a proprietary cleaning agent and follow the manufacturer's instructions. The heating/cooling block is made of aluminium, therefore an agent such as Neutracon (from Decon Labs Ltd.), suitable for nonferrous metals should be used. However, remember that other parts of the unit are made of ferrous materials and may be damaged by spillage onto them.

HEATER BLOCK REMOVAL

 Removing the block may expose the possibility of electric shock; disconnect the unit from the mains and allow to cool below 50°C before continuing. Remove the four screws on the underside of the unit and drop the block/fan assembly down. Unplug the assembly from the wiring and remove the assembly from the unit. Refitting the block/fan assembly is the reverse procedure; be careful of the wires as you refit the block.

*Underside of the TC-3000
(TC-3000X is identical)*

Four screws which hold the
block/fan assembly in place

CLEANING YOUR CASE

The outer case of the TC-3000 and TC-3000X can be cleaned with a cloth dipped in water or ethanol (50:50 methanol/water or 99:1 water/formaldehyde can also be used). No part of the case or cover should be immersed in the solvents. Do not use aggressive solvents such as acetone, or abrasive cleaners. Before using any cleaning or decontamination method except those recommended here, the responsible person should check with Bibby Scientific Ltd that the proposed method will not damage the equipment.

SOME FREQUENTLY ASKED QUESTIONS

Q1 What is the 'Pause' function at the start of the program used for?

- A1** Some users prefer to preheat the heated lid before placing the samples into the unit. The pause feature is used to stop the unit after the 4 minute heated lid preheat step. It will also sound an audible alarm indicating that the machine is ready for the sample tubes or plate to be added. Pressing the "Pause" key followed by the "Enter" key will commence the remaining program.

Q2 Why do you have to press some keys twice to make the unit do what you want it to do?

- A2** With many of the program features there is a fail-safe function which asks if you are sure you want to perform a particular action. You are required to press one key to select a function followed by a second to accept it.

Q3 What is 'Hot Start'?

- A3** A "hot start" is used for reducing non-specific product formation during the first cycle of amplification. The 'Hot Start' programming step is used to pause the instrument at a specific temperature, typically around 70°C, after the initial template denaturation. The reason is to allow the manual addition of unmodified Taq DNA polymerase which may lose activity if added during the initial 5min denaturation. Heat-activated Taq or Hot Start enzymes do not require this step.

Q4 What is the incremental function?

- A4** Incremental timing and temperature are used to increase or decrease either the time or temperature incrementally over the number of cycles in a stage.

Incrementation of extension time is used with 'Long PCR' which is when large template fragments are to be amplified (e.g. 27kb lambda DNA, 40kb genomic DNA).

Decremental temperature is used for protocols such as 'Touchdown PCR' where one starts with a high annealing temperature in the first cycle and gradually decreases the temperature over the number of cycles in the stage. This ensures that only the specific product is amplified.

Q5 The heat-up ramp-rate in the specification is 3.6°C/second yet I can only manually set it to 2.0°C/second. Why is this?

- A5** If you set the ramp-rate to max the block will heat at 3.6°C/second the unit sets its own control parameters. If you program a ramp-rate the control parameters reduce the ramp-rate. This is a feature of the software and not the actual ramp-rate.

If you require further technical or applicational assistance please contact Bibby Scientific Ltd at: **technehelp@bibby-scientific.com**.

For servicing information please contact: **service@bibby-scientific.com**.

We are continually striving to improve our thermal cyclers and software. If you have any comments and suggestions on how we can do things better please send them to us at:

techne@bibby-scientific.com

ADDITIONAL INFORMATION

Brief fault finding notes and a list of replacement parts are given in this section.

Note that this equipment should only be dismantled by properly trained personnel. Removing the outer cover exposes potentially lethal mains voltages. There are no user serviceable parts within this equipment.

FAULT FINDING

Should you have any problems with your unit which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Please include details of the fault observed and remember to return the unit in its original packing. Bibby Scientific Ltd accepts no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier, giving the full serial number of the unit and software version number (shown when the unit is first switched on).

FUSES

If neither the power light nor display on the front panel is lit, one of the two fuses may have blown. Check that there is no external cause, such as a faulty plug or lead. Check both fuses and replace the faulty fuse with a new one of the correct value (fuse values are given on the label next to the power inlet). Note that fuses should only be replaced by a qualified electrician.

The holder for the two fuses is built into the mains input socket. First remove the power cable and then gently prise the fuse drawer open with a flat-bladed screwdriver or similar tool.

Each fuse can be removed by using the screwdriver as a lever.

Exchange both fuses in the fuse holder for a working fuse of the correct value. Finally, replace the fuse drawer in the fuse compartment and push the drawer shut.

Fuses which blow repeatedly indicate a serious fault and you should return the unit to your supplier for repair.

THE HEATED LID OVER-TEMPERATURE CUT-OUT

The heated lid is fitted with an independent circuit to protect it from overheating. In the unlikely event of an over-temperature problem with the lid, the unit is fitted with a thermal fuse which will remove power to the heater plate should the maximum temperature be exceeded. If this fault occurs, contact your local supplier for repair.

INSULATION TESTING

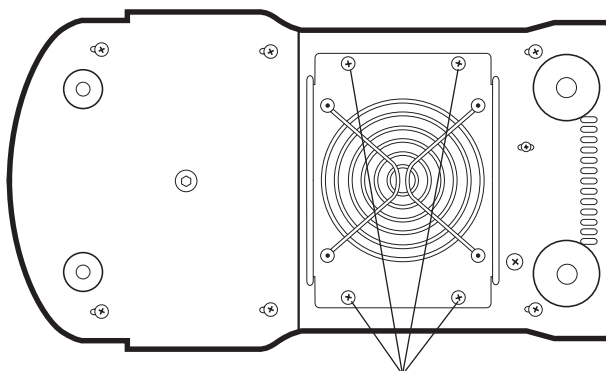
This equipment is fitted with RFI suppression circuitry. Any check of the electrical insulation by means of high voltage dielectric testing (for example as in BS EN 61010-1) must be carried out using only a DC voltage.

This unit contains semiconductor components which may be damaged by electric field effects.

INTERCHANGEABLE BLOCKS

The block can be removed and replaced by another of the same sort or of a different sort. The software will register the change and set up the unit for the new block.

Remove the four screws on the underside of the unit and drop the block/fan assembly down. Unplug the assembly from the wiring and remove the assembly from the unit. Refitting the block/fan assembly is the reverse procedure; be careful of the wires as you refit the block.

Underside of the TC-3000
(TC-3000X is identical)

Four screws which hold the
block/fan assembly in place

ACCESSORIES

The following accessories can be obtained from Bibby Scientific Ltd or your Bibby Scientific Ltd dealer:

Item No	Description	Quantity
FTC3/02/B	0.2ml Tube Block for TC-3000	1
FTC3/05/B	0.5ml Tube Block for TC-3000	1
FTUB02TW	0.2ml Micro tubes	Pack of 1000
FTUB05TW	0.5ml Micro tubes thin wall	Pack of 1000
FTC3X/02/B	0.2ml Tube block for TC-3000X	1
FTC3X/05/B	0.5ml Tube block for TC-3000X	1
FSTRIP2	Strips of 0.2ml tubes for use with flat or domed caps	Pack of 125 strips
FSTRIP3	Strips of 0.2ml tubes, each tube has individual flat cap	Pack of 120 strips
FSTRIP4	Strips of 0.2ml tubes with attached, hinged strip of domed caps	Pack of 125 strips
F8CAP2	Strips of 8 domed caps for use with FSTRIP2	Pack of 125 strips
F8CAP3	Strips of 8 flat caps for use with FSTRIP2	Pack of 125 strips
FCOOL	Mini cooler. Suitable for all 0.2ml tubes, strip tubes & plates	1
F24PLNS	24 well plate, non-skirted	40 plates
F48PLNS	48 well plate, non-skirted	20 plates
F24MAT	24 well sealing mat (silicone) for use with F24PLNS	10 mats
F48MAT	48 well sealing mat (silicone) for use with F48PLNS	10 mats

REPLACEMENT PARTS

The following replacement parts can be obtained from Bibby Scientific Ltd or your Bibby Scientific Ltd dealer:

Item No	Description	Quantity required
FCABRTUK	Mains cable and plug, 230V UK	1
FCABRTEU	Mains cable and Schuko plug, 230V Europe	1
FCABRTUS	Mains cable and 3-pin plug, 100V/120V US type	1
6501231	Fuse T2.5A (TC-3000 230V units)	2
6500129	Fuse T4A (TC-3000 120/100V units)	2
6500342	Fuse T6.3A (TC-3000X)	2
6106196	Heated lid mat (TC-3000X)	1

**Bibby Scientific Ltd**

Beacon Road · Stone
Staffordshire ST15 0SA
United Kingdom
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 813748
email: techne@bibby-scientific.com
www.bibby-scientific.com

Bibby Scientific France SAS

ZI du Rocher Vert - BP 79
77793 Nemours Cedex
France
Tel: +33 1 64 45 13 13
Fax: +33 1 64 45 13 00
e-mail: bsf@bibby-scientific.fr
www.bibby-scientific.com

Techne Inc.

3 Terri Lane Suite 10
Burlington NJ 08016
USA
Tel: 800-225-9243
Fax: 609-589-2571
email: labproducts@techneusa.com
www.techneusa.com

Bibby Scientific Italia Srl

Via Alcide de Gasperi 56
20077 Riozzo di Cerro al Lambro
Milano Italia
Tel: +39 (0)2 98230679
Fax: +39 (0)2 98230211
e-mail: marketing@bibby-scientific.it
www.bibby-scientific.it

Bibby Scientific (Asia) Ltd

Room 607 Yen Sheng Centre
64 Hoi Yuen Road Kwun Tong
Kowloon Hong Kong
Tel: +852 3583 1581
Fax: +852 3583 1580
e-mail bibby@bibby-scientificasia.com
www.bibby-scientificasia.com