

# BD Accuri<sup>™</sup> C6 Software User Guide

Science is hard. Flow cytometry should be easy.  $^{^{\text{TM}}}$ 



#### **TABLE OF CONTENTS**

1		ODUCTION TO BD ACCURI C6 SOFTWARE	
	1.1	Starting BD Accuri C6 Software	
	1.2	BD Accuri C6 Software Workspace	
	1.3	Opening a New BD Accuri C6 Software Workspace	
	1.4	Exiting BD Accuri C6 Software	
	1.5	Using the Example BD Accuri C6 Software File	3
2	VALI	DATING THE PERFORMANCE OF THE BD ACCURI C6 FLOW CYTOMETER	5
	2.1	Running Validation Beads (Daily Start-Up)	
	2.2	Saving Validation Bead Data	
	2.3	Analyzing and Recording Validation Bead Data	
	2.4	Monitoring Validation Bead Data	
	2.5	Troubleshooting Validation	
2		ACQUISITION	16
3			
	3.1 3.2	Viewing the <i>Collect</i> Tab	
	3.2	Collecting Sample Data	
		3.2.1 Setting the Fluidics Rate	
		3.2.2 Setting the Threshold	
		3	
		3.2.4 Naming the Sample	
		3.2.5 Running the Sample	
		3.2.7 Pausing Data Collection	
	3.3	Ending a Data Collection Session	
	3.4	Creating Plots	
	3.5	Changing Plot Specifications	
	3.6	Using Gates and Markers	
	3.0	3.6.1 Creating a New Gate	
		3.6.2 Applying a Gate to a Plot	
		3.6.3 Creating and Applying Nested Gates	
	3.7	Moving and Resizing Regions	
	3.8	Changing the Number of Events in a Plot	
	3.9	Naming Plot Axes	
	3.10	Zooming on a Plot	
	5.10	3.10.1 Basic Zoom	
		3.10.2 Zooming to a Specified Channel Range	
	3.11	Saving a BD Accuri C6 Software File	
	0.11	3.11.1 Auto-Saving Files	
		3.11.2 Manually Saving Files	
	3.12	Creating a BD Accuri C6 Software Template	43
		Understanding Fluorescence Spillover	
	0.10	3.13.1 Recognizing Fluorescence Spillover	
		3.13.2 Correcting Fluorescence Spillover	
		3.13.3 Troubleshooting Color Compensation	
	3.14	Changing Parameters	
	3.15	Copying and Pasting Plots	
	3.16	Printing Data	
	3.17	Exporting and Importing Files	
4	ANAL	YZING SAMPLE DATA	54

	4.1 4.2		ng the <i>Analyze</i> Tabg up Plots	
	4.2		Copying Plots from the <i>Collect</i> Tab	
		4.2.2	Creating Plots	57
			Creating an Overlay Histogram	
	4.3	Viewir	ng Plots	59
5	CREA	TING A	A STATISTICS TABLE	61
	5.1		ng the Statistics Tab	
	5.2		ng the Master Statistics Table	
	5.3		ewing a Plot in the Statistics Tab	
	5.4	Copyi	ng Data into Other Applications	64
6	MAIN		IG THE BD ACCURI C6 FLOW CYTOMETER	
	6.1		ing the SIP	
	6.2		ing the Flow Cell	
			Running the Unclog Cycle	
	6.2		Running an Extended Clean of the Flow Cell	
	6.3 6.4		ing the Fluidics Linestaminating the Fluidics System	
	6.5		the BD Accuri C6 Flow Cytometer for Precise Volume Measurements	
_		_	-	
7			PEAK POSITION WITH VIRTUALGAIN™	
	7.1 7.2		ng VirtualGain ng VirtualGain	
	7.2		ving VirtualGainving VirtualGain	
	PENDI		BD ACCURI C6 SOFTWARE MENU QUICK REFERENCE	
API	PENDI	хв	ADVANCED FLUIDICS SETTINGS	78
APF	PENDI	_	TRACKING USER ACTIVITY	
	C.1	Install	ing the User Tracking Module	80
	C.2		the Tracking Feature for the First Time	
	C.3		g, Deleting, and Modifying User Accounts	
			Adding User Accounts  Deleting User Accounts	
			Changing a Password	
	C.4		g In and Signing Out	
	C.5		ring a Forgotten Administrator Password	
	C.6		oring User Activity	
API	PENDI	X D	BD ACCURI C6 ANALYSIS SOFTWARE	85
۸Ы	PENDI	v =	SELECTABLE LASERS	96
AFI	E.1		ing the Selectable Lasers Module	
	E.2		ting Proper Function After Installation	
	E.3		ating Selected Laser Configuration	
	E.4	Optica	al Filter Placement	92
	E.5		able Laser Application Examples	
			2 Blue 2 Red Configuration Examples	
		E.5.2	4 Blue Configuration Examples	97
APF	PENDI		NHANCED ANALYSIS FEATURES	
	F.1		ng a Live Gate	
	F.2	Renar	ning Plots and Regions	101

F.3	Colori	ing Events in a Region	102
		ing Publication-Quality Images of Plots	
		zing Batches of Samples	
	F.5.1	Viewing the Batch Analysis Tab	103
	F.5.2	Running a Batch Analysis	104
		Exporting Data	
APPENDI	IX G	FCS KEYWORDS	106
APPENDI	IX H	EXAMPLE .CSV FILE	109
APPENDI	IX I T	FECHNICAL SUPPORT	110
For T	echnica	al Support, contact	110

7820095-01 Rev-0 iii

#### **TABLE OF FIGURES**

Figure 1-1. Collect Tab Workspace	2
Figure 1-2. New BD Accuri C6 Software Workspace	3
Table 1-1. Experimental Design for HPB 4 Color Tutorial File	
Figure 2-1. Open the Bead Template	
Figure 2-2. Select an Empty Well	
Figure 2-3. Run Limits: Disable Run Unlimited	
Figure 2-4. Run Limits: 50000 Events	
Figure 2-5. Sample Name: 8-Peak Beads	
Figure 2-6. Select an Empty Well	
Figure 2-7. Run Limits: 50000 Events	
Figure 2-8. Sample Name: 6-Peak Beads	
Figure 2-9. Run Limits: 2 Minutes	
Figure 2-10. Plot with Bead Doublets	
Figure 2-11. Gate Applied to 8-Peak Bead Plot	10
Figure 2-12. Zoomed View of Plot	
Figure 2-13. Percentage of a population within a given marker or region appears on plots; the	
percent CV value appears only in Statistics tables of BD Accuri C6 Software	
Figure 2-14. 6-Peak Validation Bead Data	
Figure 2-15. Statistics Tab: 8-Peak Bead Data from Successive Days	
Figure 3-1. Collect Tab Workspace	
Table 3-1. Collect Tab Controls	
Figure 3-2. New BD Accuri C6 Software Workspace	
Table 3-2. Suggested Starting FSC-H Threshold Settings for Various Cell Types	
Figure 3-3. Threshold Settings Warning Message	21
Figure 3-4. Primary Threshold Drop-Down List	
Figure 3-5. Threshold Settings Dialog Box	
Figure 3-6. Run Limits Controls	
Figure 3-7. Sample Name Field	
Figure 3-8. BD Accuri C6 Software Workspace after Collecting a Sample	
Figure 3-9. BD Accuri C6 Software Workspace with Empty Sample Well and Empty Plots	
Figure 3-10. View of a Plot	
Figure 3-11. New Density and Histogram Plots	
Figure 3-12. Set Plot Specs Dialog Box	
Figure 3-13. Using Polygonal Gating Tool	
Figure 3-14. Using the Vertical Marker	
Figure 3-15. Using the Horizontal Marker	
Figure 3-16. Selecting a Gating Option	
Figure 3-17. Plot Gated to Include P1	
Figure 3-18. First Gate for Creating Nested Gates, Drawn in Plot 4	
Figure 3-19. Applying the Parent Gate to Plot 5	
Figure 3-20. Parent Gate Applied to Dot Plot	
Figure 3-21. Second Gate for Creating Nested Gates	
Figure 3-22. Applying the Child Gate	
Figure 3-23. Third Plot with Nested Gate Applied (R1 in P1)	36
Figure 3-24. Statistics of Plot with Nested Gate	
Figure 3-25. Region P1 in the plot has been selected, as indicated by bolded outline	
Figure 3-26. Moved region in the <i>Batch Analysis</i> tab	
Figure 3-27. Before and After Changing Events Displayed	
Figure 3-28. Events Display Settings Dialog Box	
Figure 3-29. Plot with Events Display Settings Applied	

Figure 3-30. X-Axis Label	39
Figure 3-31. Rename Parameters Dialog Box	39
Figure 3-32. Rename Parameters Dialog Box with Axis Label Drop-Down List	40
Figure 3-33. Before and After Using Zoom Tool	40
Figure 3-34. Plot Spec Dialog Box: Set Min and Max Channel Values for the X- and Y-Axes.	41
Figure 3-35. Title Bar with File Name	42
Figure 3-36. Auto-Save Settings Dialog Box	
Figure 3-37. Save BD Accuri C6 Software File	
Figure 3-38. Save BD Accuri C6 Software Template	
Figure 3-39. Fluorescence Spillover in Different Plots	
Figure 3-40. Corrected Fluorescence Spillover	
Figure 3-41. Placing a Quadrant Tool	
Figure 3-42. Statistics Table Displaying Median Values	
Figure 3-43. Compensation Settings Dialog Box	
Table 3-3. Example of Fluorochrome Spillover per Channel,	
Figure 3-44. Results of Subtracting Spillover	
Figure 3-45. Plot Displaying Overcompensation	
Figure 3-46. Set Plot Specs Dialog Box with Hide 1st decade Disabled	
Figure 3-47. Properly Compensated Data after First Decade Is Rehidden	
Figure 3-48. Set Plot Drag and Drop Format Dialog Box	
Figure 3-49. Plots Selected for Printing	
Figure 3-50. Open an FCS file	
Figure 4-1. Analyze Tab Workspace	
Table 4-1. Analyze Tab Controls	
Figure 4-2. Selecting Plots to Copy from the Collect Tab	
Figure 4-3. Plot List Containing Copied Plots	
Figure 4-4. Blank Overlay Histogram Plot	
Figure 4-5. Overlay Histogram Plot with Data	
Figure 4-6. Overlay Histogram Legend	
Figure 4-7. Overlay Histogram Legend with Color Palette	
Figure 4-8. Analyze Tab: Plots with Gating Applied Figure 4-9. Analyze Tab: Two Samples Displayed with the Same Plots	
· · · · · · · · · · · · · · · · · · ·	60 61
Figure 5-1. Statistics Tab Workspace	61 62
Figure 5-1. Statistics Tab Controls	62 62
Figure 5-3. Creating Master Statistics Table: Adding Prots	62 62
Figure 5-4. Plot Preview	
Figure 5-5. Sample Selector List and Plot	บว
Figure 7-1. Before and After Applying VirtualGain	
Figure 7-2. Setting Up Histograms for VirtualGain	03 70
Figure 7-3. Aligning Plots	 71
Figure 7-4. Pick the Sample to Align	 71
Figure 7-5. Move the Peak Definition Marker	 72
Figure 7-6. Align Additional Samples	
Figure 7-7. Black Asterisk Identifier	
Figure 7-8. Plot with Black Asterisk	
Figure 7-9. VirtualGain Applied in an Overlay Histogram	
Figure 7-10. Toggle between VirtualGain Applied (Left) and Not Applied (Right)	
Table A-1. BD Accuri C6 Software Menus	
Figure B-1. Custom Radio Button	
Figure B-2. Set Core Size	

7820095-01 Rev-0 v

Table B-1. Core Size/Flow Rate Combinations	79
Figure C-1. User Tracking Installer Icon	80
Figure C-2. Installation Dialog Box	80
Figure C-3. Username and Password Dialog Box	81
Figure C-4. The Users Information Box	81
Figure C-5. Adding a New User	82
Figure C-6. Deleting a User	82
Figure C-7. Username and Password Dialog Box	83
Figure C-8. userUsage Log	84
Figure C-9. CytometerSupportFiles Folder	84
Figure E-1. Selectable Lasers Installer	86
Figure E-2. Install Wizard: Choose Activation Keys Directory	87
Figure E-3. Install Wizard: Install the Module	88
Figure E-4. Install Wizard: Setup Completed Successfully	88
Figure E-5. Selectable Laser Controls in the BD Accuri C6 Software Workspace	
Figure E-6. Selectable Laser Options: 3 Blue 1 Red	89
Figure E-7. Selectable Laser Options: 2 Blue 2 Red	90
Figure E-8. Selectable Laser Options: 4 Blue	90
Figure E-9. Evaluation of 6-Peak Validation Beads for the 2 blue 2 red Configuration	
Figure E-10. Evaluation of 8-Peak Validation Beads for the 4 Blue Configuration	
Figure E-11. Renamed Data Well to Indicate 4 Blue Configuration	
Table E-1. 3 Blue 1 Red:    Configuration 1 (Standard Filters)	
Table E-2. 3 Blue 1 Red: Configuration 2	
Table E-3. 2 Blue 2 Red Configuration	
Table E-4. 4 Blue: Configuration 1	
Table E-5. 4 Blue: Configuration 2	
Table E-6. 4 Blue: Configuration 3	
Table E-7. Detector and Filter Configuration for 2 Blue 2 Red, Examples 1 and 2	
Figure E-12. Gating Example Using the 2 Blue 2 Red Configuration	
Figure E-13. BD Accuri C6 Software (PN 653126).	
Table E-8. Detector and Filter Configuration for Selectable Lasers 4 Blue Example 1	
Figure E-14. 4 Blue Configuration and the Optional 780/60 Optical Filter at Detector FL3	
<b>Table E-9.</b> Detector and Filter Configuration for Selectable Lasers 4 Blue, Example 2	
Figure E-15. HPB Stained with CD45-FITC, CD4-PE, CD8-PE-Texas Red and CD3-PE-Cy5.	
Figure F-1. Creating a Live Gate	
Figure F-2. Enable Renaming of the Plot	
Figure F-3. Type a New Plot Name	
Figure F-4. New Plot Name	
Figure F-5. Select a Region to Color	
Figure F-6. Select a Color for a Region	
Figure F-7. Colored Events with a Region	
Figure F-8. Batch Analysis Tab Workspace	
Table F-9. Batch Analysis Tab Controls	
Figure F-10. Export Buttons in the Batch Analysis Tab	
Table G-1. Standard FCS Tags	
Table G-2. Custom FCS Tags	
Table G-3. FCS Tags Not in BD Accuri C6 Software Exported Files	ıUÖ

#### 1 INTRODUCTION TO BD ACCURI C6 SOFTWARE

BD Accuri C6 Software controls the BD Accuri C6 flow cytometer system in order to acquire data, generate statistics, and analyze results.

BD Accuri C6 Software provides the following features:

- Tabbed views for collection, analysis, and statistics
- Digital signal processing and color compensation at any time
- Drag and drop plots
- Export of files in FCS 3.0 format
- Seamless data file importation into FCS Express
- Enhanced Analysis upgrade adds:
  - Drag and drop of publication-quality images, event coloring, live gating
  - Batch analysis of sample data

#### 1.1 Starting BD Accuri C6 Software

To open BD Accuri C6 Software:

• Double-click on the *Accuri C6 Software* icon on the computer desktop. BD Accuri C6 Software opens a new, blank workspace.

**NOTE:** If BD Accuri C6 Software displays the message *Extra startup time needed* due to cleaning or improper shutdown, the cytometer will take several more minutes than usual to recover and return to the green-light ready state. This may occur on initial cytometer startup after installation. It will also occur after an interruption of power to the unit.

#### 1.2 BD Accuri C6 Software Workspace

The main BD Accuri C6 Software window is called the BD Accuri C6 Software workspace. The workspace contains controls and displays that provide access to all functions required for data acquisition and analysis. The workspace is organized on four separate tabs:

- Collect—Contains controls for setting up data collection and acquiring data (see chapter 3 for details).
- Analyze—Allows analysis of multiple samples simultaneously (see chapter 4 for details).
- Statistics—Displays statistical information (see chapter 5 for details).
- Batch Analysis—Contains controls for analyzing batches of sample data and is available with the Enhanced Analysis upgrade (see APPENDIX F for details).

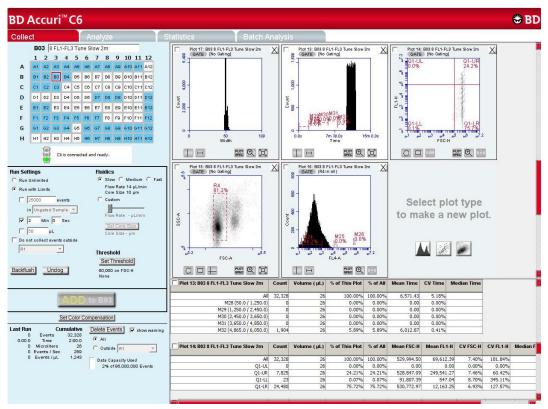


Figure 1-1. Collect Tab Workspace

## 1.3 Opening a New BD Accuri C6 Software Workspace

A new workspace (Figure 1-2) displays a single FSC-A versus SSC-A density plot that is pre-zoomed to channel values of 1,600,000 and 800,000 respectively. Run settings will be set to *Run Unlimited*, the threshold will be set to channel 80,000 on the FSC-H signal, and no color compensation values are set.

A new workspace can be used to create an analysis template and to collect a new dataset.

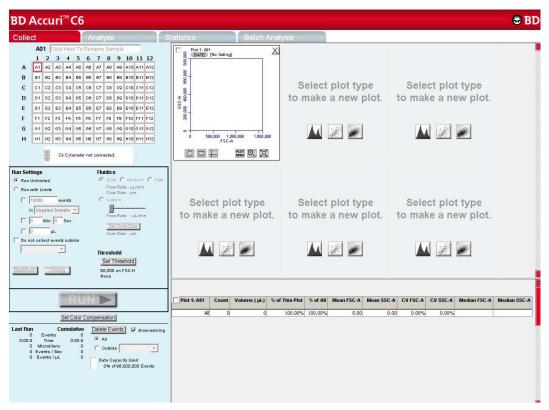


Figure 1-2. New BD Accuri C6 Software Workspace

To open a new workspace:

Do one of the following:
 If BD Accuri C6 Software is not already open, double-click on the BD Accuri C6 Software icon on the computer desktop. If BD Accuri C6 Software is already open, select File > New Workspace. If desired, save any unsaved changes to the previous workspace when prompted.

# 1.4 Exiting BD Accuri C6 Software

- Select File > Quit.
- If prompted to save changes to the BD Accuri C6 Software workspace, do one of the following:
  - a. Click on the Yes button to save changes.
  - b. Click on the *No* button to close the software without saving changes.
  - c. Click on the Cancel button to cancel the exit and keep the software open.

# 1.5 Using the Example BD Accuri C6 Software File

An example BD Accuri C6 Software file of a four-color analysis of human peripheral blood (HPB 4 Color Example.c6) can be downloaded from the BD Accuri website (<a href="www.accuricytometers.com/resources/tutorials/">www.accuricytometers.com/resources/tutorials/</a>). This data file can be used to explore various tools in BD Accuri C6 Software without the worry of corruption or loss of experimental data. Figures throughout this user guide show data from the example file.

To create the example file, four sample tubes were used to assess the CD3<sup>+</sup>CD4<sup>+</sup> and CD3<sup>+</sup>CD8<sup>+</sup> cell populations. These samples were prepared by staining peripheral blood with directly conjugated antibodies, followed by red cell lysis, according to standard methods.

The following table describes the experimental design:

- Tube 1—Background control (unstained)
- Tube 2—White blood cell gating control (CD45<sup>+</sup>)
- Tube 3—T-cell gating control
- Tube 4—Experimental sample

Table 1-1. Experimental Design for HPB 4 Color Tutorial File

Fluorochrome	Sample Tube			
	Tube 1	Tube 2	Tube 3	Tube 4
FITC	unstained	unstained	CD3	CD3
PE	unstained	unstained	unstained	CD4
PE-Cy7	unstained	CD45	CD45	CD45
APC	unstained	unstained	unstained	CD8

This experimental design does not contain all single-stained fluorescence controls, but only those required for determining color compensation settings to correct fluorescence spillover. For details on correcting fluorescence spillover, see section 3.13.

# 2 VALIDATING THE PERFORMANCE OF THE BD ACCURI C6 FLOW CYTOMETER

Perform a validation of the system at least once each day of use. This ensures that the cytometer is working properly before running experimental samples.

Using the same BD Accuri C6 Software file each day allows identification of trends over time. When collecting validation data, advance to the next empty well in row A-D (for 8-peak beads) or E-H (for 6-peak beads). Include the date in the Sample Naming Field for each day of validation. Start a new validation BD Accuri C6 Software bead file when all data wells are full.

#### Reagents required:

- Spherotech 8-Peak Validation Beads (PN 653144, supplied with the initial BD Accuri C6 flow cytometer shipment)
- Spherotech 6-Peak Validation Beads (PN 653145, supplied with the initial BD Accuri C6 flow cytometer shipment)
- Sheath fluid: Filtered (0.2 µm filter), deionized water plus Bacteriostatic Concentration Solution (PN 653156, supplied with the initial BD Accuri C6 flow cytometer shipment)

# 2.1 Running Validation Beads (Daily Start-Up)

#### Setup:

Verify that the BD Accuri C6 Software 8 and 6 Peak Bead Template has been copied to the computer workstation. The file is located on the BD Accuri C6 Software CD or flash drive and is on the BD Accuri website (<a href="https://www.accuricytometers.com/resources/templates/">www.accuricytometers.com/resources/templates/</a>).

Open BD Accuri C6 Software.

Select File > Open Workspace or Template.

In the Open dialog box, browse to the location of the bead template file and open the file. If adding to an established BD Accuri C6 Software bead file, browse to the location of the file.

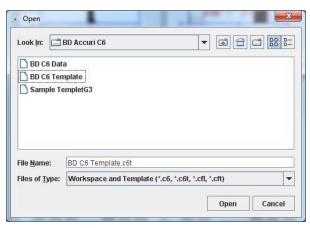


Figure 2-1. Open the Bead Template

Click on the first empty well in rows A-D to advance to the well.

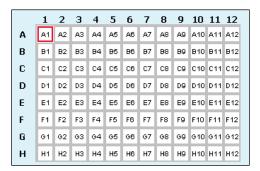


Figure 2-2. Select an Empty Well

Place a tube with 2 mL of filtered, deionized water on the SIP.

Enable the Run with Limits radio button in the Instrument Control Panel.

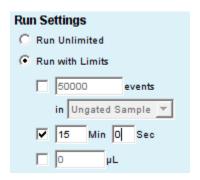


Figure 2-3. Run Limits: Disable Run Unlimited

Enable the *Time* check box next to the Min and Sec fields in the Instrument Control Panel and type in a run time of fifteen minutes.

Select the Fast radio button in the Fluidics section of the Control Panel.

Click on the RUN button to rinse the SIP.

When prompted, save the file.

Once the run is finished, click on the *Delete Sample Data* button to delete data collected during the rinse.

#### Run 8-Peak Validation Beads:

Disable the *Time* check box next to the *Min* and *Sec* fields, enable the *Events* check box and enter *50,000* into the *Events* field

Select *Ungated Sample* from the associated drop-down list.

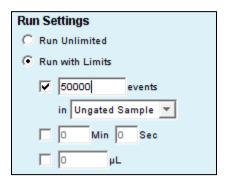


Figure 2-4. Run Limits: 50000 Events

Vortex a sample tube containing suspended 8-peak validation beads, prepared according to the package instructions. Place the tube on the SIP.

Select the Slow radio button in the Fluidics section of the Control Panel.

Click on the *RUN* button to start acquisition. Acquisition automatically stops after 50,000 total events are acquired.

**CAUTION:** Make sure the well in BD Accuri C6 Software is empty before starting the run. If the button displays *ADD TO*, the well already contains data.

**NOTE:** The R1 region may not encompass the main population of bead events on the FSC-H vs. SSC-H plot. This is common and acceptable at this stage.

Name the sample by typing a name in the text box just above the Sample Grid. Include the date in the sample name to differentiate it from samples collected on other dates.

**NOTE:** Samples can be named before, during, or after collection.

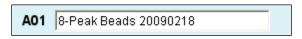


Figure 2-5. Sample Name: 8-Peak Beads

When the collection is finished, remove the sample tube and wipe off the end of the SIP with a lint-free tissue (or similar material) to minimize sample carryover.

#### Run 6-Peak Validation Beads:

Vortex a tube of suspended 6-peak validation beads, prepared according to the package instructions. Place the tube on the SIP.

Click on the first empty well in rows E-H to advance to the well.

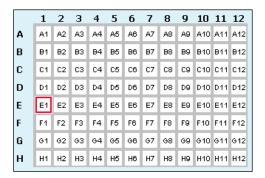


Figure 2-6. Select an Empty Well

Verify that the check box by *Events* is still enabled and set at 50000 and that *Ungated Sample* is still selected from the drop-down list.

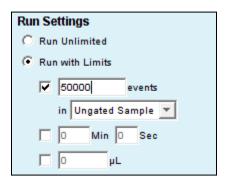


Figure 2-7. Run Limits: 50000 Events

Click on the RUN button.

**NOTE:** The R2 region may not encompass the main population of bead events on the FSC-H vs. SSC-H plot. This is common and acceptable at this stage.

Name the sample with a name that includes the date processed.



Figure 2-8. Sample Name: 6-Peak Beads

#### **End The Procedure:**

When the collection is finished, remove the sample tube from the SIP and wipe off the end of the SIP with a lint-free tissue.

Place a tube with 2 mL of filtered, deionized water on the SIP and advance to any empty data well.

Select the *Time* check box (*Min Sec*) in the Instrument Control Panel and set it for two minutes.

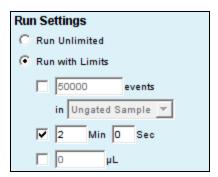


Figure 2-9. Run Limits: 2 Minutes

Click on the RUN button.

When the run is finished, leave the tube on the SIP.

#### 2.2 Saving Validation Bead Data

By default, BD Accuri C6 Software automatically saves data at the end of each sample run. Data can also be saved manually at any time. For information on saving data, see section 3.11.

#### 2.3 Analyzing and Recording Validation Bead Data

After bead data is collected, analyze the data using the *Collect* tab of BD Accuri C6 Software to ensure that the cytometer is functioning properly:

Click on the well that contains the most recent 8-peak bead data (in rows A-D).

On the first FSC-H vs. SSC-H plot (scatter plot) in the bead file, adjust the pre-drawn region (R1) to encompass the main population by dragging the border of the region (see Figure 2-10). R1 should contain at least 80% of all events.

**NOTE:** There is usually a "shadow" population (called *bead doublets* or *clumps*) that is slightly higher in FSC-H than the main cluster of beads; this is normal for these beads. Do not include the shadow group in R1.

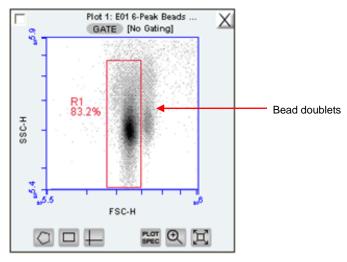


Figure 2-10. Plot with Bead Doublets

Verify that the next three plots (FL1-H, FL2-H, and FL3-H) are gated on scatter region R1 and that the plots display the message *R1 in all* next to the *GATE* button (Figure 2-11). If it is not displayed, click on the *GATE* button and select *R1 on all* events from the pop-up dialog box. For information about gates, see section 3.6.

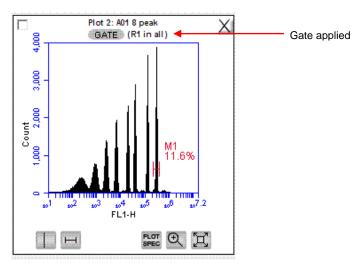


Figure 2-11. Gate Applied to 8-Peak Bead Plot

 Measure the CV of the top (brightest, far right) peak on each of the three fluorescence plots. To place the pre-drawn horizontal marker tightly around the peaks:

Use the Zoom Tool \( \bigsize \) in the plot to zoom in on the top peak (see section 3.10). Adjust the marker tightly around the peak by clicking on the marker and dragging its edges.

Click on the Expand Tool I to zoom back out.

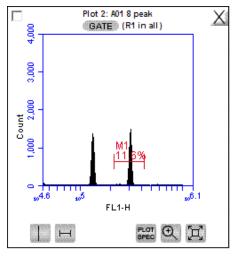


Figure 2-12. Zoomed View of Plot

Compare the bead run to the manufacturing results sent with the cytometer. If the cytometer is performing properly, the data plots should look similar to the 8-peak bead plots in Figure 2-13. Look for the following:


One main population of beads on FSC-H vs. SSC-H (a shadow population is acceptable)

Eight discernible peaks on FL1-H and FL2-H

At least six peaks on FL3-H

Top Peak CV's are less than 5%.

**NOTE:** FL4-H performance is validated with the 6-peak beads in steps 6-9.

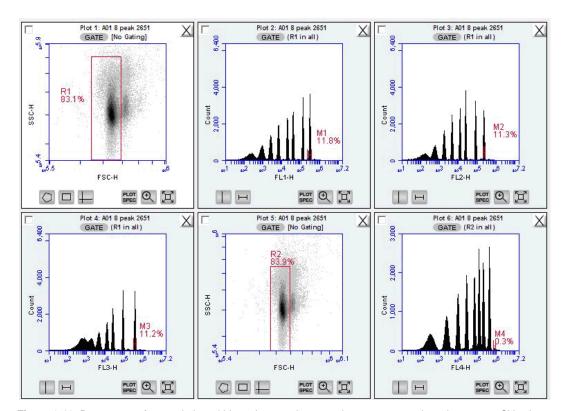

**Figure 2-13.** Percentage of a population within a given marker or region appears on plots; the percent CV value appears only in Statistics tables of BD Accuri C6 Software.

Select the well containing the most recent 6-peak bead data (in rows E-H).

Adjust the pre-drawn region R2 in the 6-peak scatter plot FSC-H vs. SSC-H (plot 5) to encompass the main population (similar to the procedure for the 8-peak beads). This population should look like an exclamation point. The R2 region should encompass the entire exclamation point (Figure 2-14).

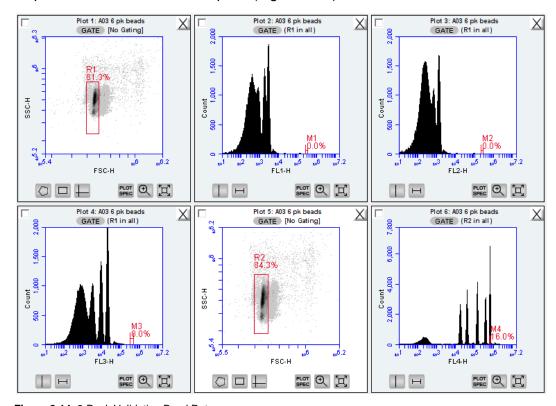


Figure 2-14. 6-Peak Validation Bead Data

Verify that the FL4-H plot is gated on region R2. If it is not, click on the *GATE* button and select *R2 on all events* from the pop-up dialog box.

• Measure the CV of the top (brightest, far right) peak by adjusting the marker in the FL4-H plot so that it is placed tightly around the peak. See the plots in Figure 2-14 for an example of acceptable 6-peak bead data. Look for the following:

One main population of beads on FSC-H vs. SSC-H

Six peaks on FL4-H

If desired, record the following information for each parameter in the BD Accuri C6 Flow Cytometer Log (available on the BD Accuri C6 Software CD or flash drive):

Number of peaks

Mean channel numbers for the top peaks and forward scatter

CVs for the top peaks and forward scatter

If this is the first validation run, send a copy of the file to BD Accuri Technical Support for analysis. See Appendix I for email address.

### 2.4 Monitoring Validation Bead Data

If a single BD Accuri C6 Software file is used to collect validation bead data, it is easy to monitor cytometer performance over time using the *Statistics* tab.

To monitor the bead data:

- Save the 8- and 6-peak validation bead data from each day in separate wells.
   Grouping the 8-peak data runs separately from the 6-peak runs in the 96-well grid simplifies use of the Statistics tab.
- Create a table in the Statistics tab that contains the mean channel numbers and CVs for the top peaks on each detector and the forward scatter (see chapter 5 for details on creating tables in the Statistics tab).

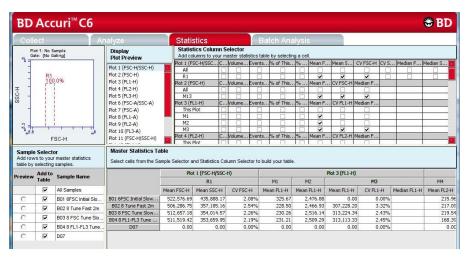


Figure 2-15. Statistics Tab: 8-Peak Bead Data from Successive Days

 Compare statistics over time for trends or sudden changes in mean values to validate the cytometer's performance. The BD Accuri C6 flow cytometer reports arithmetic means.

### 2.5 Troubleshooting Validation

The following conditions may indicate a problem with the cytometer or the validation beads:

- Very broad CV (> 5.0%) or multiple populations for FSC-H on the 8- or 6peak beads (excluding the doublet population).
- Fewer than eight peaks for FL1 or FL2.
- Fewer than six peaks for FL3 or FL4.

To troubleshoot validation:

- If bead data were not acquired using the slow rate, select *Slow*, resuspend the beads, and recollect the data.
- If the beads have been diluted for more than one week, kept at room temperature or warmer, or exposed to light for long periods of time, their performance may be substandard. Make new bead suspensions and run the bead sample again.
- There may be a bubble or clog in the flow cell. Do one or more of the following:
  - a. Run the bead sample again.

- b. Remove the tube from the SIP, place an empty tube under the SIP, and click on the *Unclog* button. When the cycle is finished (BD Accuri C6 Software displays a green Traffic Light), run the bead sample again.
- c. Remove the tube from the SIP, place an empty tube under the SIP, and click on the *Backflush* button. When the cycle is finished (BD Accuri C6 Software displays a green Traffic Light), run the bead sample again.
- d. Contact BD Accuri Technical Support (See final page of this manual for region-specific contact information.)

#### 3 DATA ACQUISITION

The *Collect* tab is used to set data collection criteria, start and stop data acquisition, and view data on collected samples. The tab contains buttons and controls for performing the following functions:

- Acquiring data
- Creating plots (histogram, dot, or density) for viewing data
- Setting stop criteria and thresholds
- Controlling the fluidics
- Using regions and markers to create gates and obtain statistics
- Printing plots and statistics
- Setting up gating strategies
- Performing fluorescence compensation
- · Importing and exporting data

#### 3.1 Viewing the *Collect* Tab

The *Collect* tab is automatically displayed when BD Accuri C6 Software opens. This tab can also be viewed by clicking on *Collect* from the *Analyze*, *Statistics*, or *Batch Analysis* tabs.

The *Collect* tab is organized into two major sections:

- Instrument Control Panel—Panel on the left side of the window that contains controls for collecting data.
- Data display—Large area on the right side of the window that shows sample data in plots and in a Statistics Table.

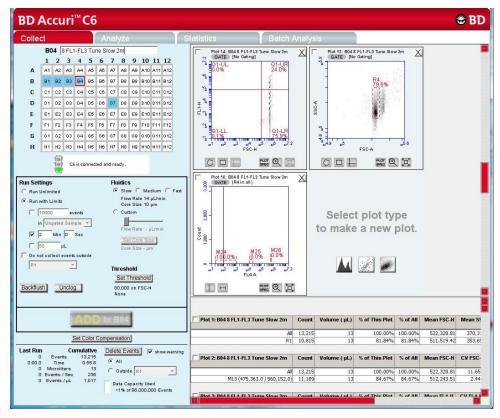


Figure 3-1. Collect Tab Workspace

The following table describes each of the controls and indicators in the *Collect* tab:

Table 3-1. Collect Tab Controls

Table 3-1. Collect Tab Collitions		
Control	Description	
Sample Naming Field	Text box for naming the current sample.	
Sample Grid	Matrix laid out in the configuration of a 96-well plate to help organize experiments and collect data from sample tubes. BD Accuri C6 Software acquires each sample into its own well in the Sample Grid. The wells can be filled with data in any order.	
	The wells are color-coded:	
	White—Does not contain data.	
	Blue—Contains data.	
	Red outline—Currently selected for viewing or collecting data.	

Control	Description
Traffic Light and Message	Indicator that displays BD Accuri C6 Software's readiness and system messages. Before data collection can begin, the software must display a green Traffic Light with the message C6 is connected and ready.
	The Traffic Light status is color-coded:
	Green—BD Accuri C6 Software is ready to collect data or is collecting data.
	Yellow—The cytometer is preparing to perform an action or a non-critical error has occurred.
	Red—A critical error has occurred.
Run Settings	Contains a set of controls that allow criteria definition for automatically stopping data collection. See section 3.2.3 for details.
Fluidics	Contains a set of controls for defining flow rate and core size. See section 3.2.1 for details.
Backflush	Performs a backflush to clear the SIP. See section 6.1 for details.
Unclog	Performs an unclog cycle to purge the flow cell of debris. See section 6.2.1 for details.
Set Threshold	Sets the event threshold to gate out debris and noise from samples. The default value is 80,000 on FSC-H. See section 3.2.2 for details on setting threshold values.
RUN/PAUSE/ADD TO	Toggle button that performs the following functions:
	RUN—Starts the sample acquisition.
	PAUSE—Pauses the acquisition. Click on ADD TO to resume data collection.
	ADD TO—Allows additional sample collection into a well that already contains data.
Set Color Compensation	Opens the Color Compensation dialog box for correcting fluorescence spillover. See section 3.13 for details.
Acquisition Counters	Displays the following information about the most recent acquisition for the selected well (Last Run) and all acquisitions for the selected well (Cumulative) in real-time:
	Events—Number of events sampled.
	Time—Elapsed acquisition time.
	Microliters—Volume of acquired sample.
	<ul> <li>Events/sec—Events acquired per second. When the run is completed, this is the average value.</li> </ul>
	• Events/µL—Events acquired per microliter. When the run is completed, this is the average value.
Delete Events	Permanently deletes all events from the current sample. Enable the <i>Show warning</i> check box to display a warning message before deleting sample data. Also contains a <i>Data Capacity Used</i> meter that displays the amount of data storage capacity currently used in BD Accuri C6 Software.  Gated events can be deleted with the <i>Enhanced Analysis Features</i> .

Control	Description
Plots Pane	Area displaying two rows of plot corrals for graphically viewing data on the selected sample. Scroll up or down to view additional plots. Each plot corral contains buttons for creating histogram, density, and dot plots. For information on creating and using plots, see section 3.4.
Statistics Table	Table below the plots that displays statistical information on individual plots.

#### 3.2 Collecting Sample Data

Events can be added to a well containing data. When a data well already contains data, the *RUN* button displays *ADD TO*. Run limits may need to be adjusted to accommodate additional data. Each data well holds a maximum of 1 million events.

BD Accuri C6 Software must display a green Traffic Light and the message *C6 is connected and ready* to collect data.

Figure 3-2 shows a new workspace. Only a density plot of linear FSC-A vs. linear SSC-A is displayed. The plot is already zoomed to show channels 0 to 1,600,000 on FSC-A and 0 to 800,000 on SSC-A.

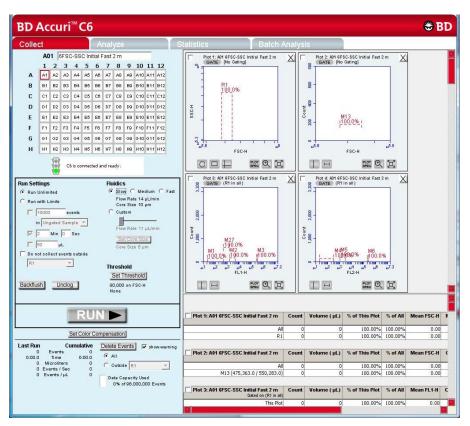


Figure 3-2. New BD Accuri C6 Software Workspace

To collect sample data:

- In the *Collect* tab, do one of the following:
  - Select a plate type and name the plate (optional).
  - · Open a template.
- Create new plots (optional, see section 3.4).
- Define acquisition settings (see sections 3.2.1 through 3.2.3).
- If necessary, acquire some events to define regions, gates, and other settings.
- Perform analysis in the Collect or Analyze tab.

#### 3.2.1 Setting the Fluidics Rate

The system can accommodate an upper limit of 10,000 events per second, but it is recommended to acquire samples at a rate of 2,500 events per second or less to ensure the best data resolution.

To set the fluidics rate:

 Click on the Slow, Medium, or Fast radio button in the Fluidics section of the Collect tab.

**NOTE:** It is recommended to start data collection on slow and observe the data rate. The setting can then be adjusted to medium or fast, if necessary. If the data rate is near or above 10,000 total events per second on the *Slow* setting, there are several possible solutions:

- Increase the primary threshold channel, taking care that the increase does not remove cells of interest from the data set.
- Include a secondary threshold, taking care not to exclude cells of interest.
- Dilute the sample.

Fluidics rate settings and sample core size can be adjusted to accommodate very small or very large particles. See APPENDIX B for more details on custom flow rates and core sizes.

### 3.2.2 Setting the Threshold

Use thresholds to gate out light scatter and/or fluorescence signals caused by debris in cell samples and electronic noise inherent in the system. When set correctly for any given sample set, resolution of particle or cell light scatter and fluorescence signals is greatly improved and data set size often can be reduced. By default, BD Accuri C6 Software is set to a primary threshold of channel 80,000 on FSC-H.

Notes on setting thresholds:

Threshold settings can be changed before, during, or after data acquisition, but
the most consistent, predictable results will be obtained if threshold settings are
chosen before *final* data collection for any given experiment. The primary
threshold is the parameter that triggers data collection. Optional secondary
thresholds can be applied to filter out additional data.

All thresholds are set on the Height signal for any given parameter. For best
results when setting or changing thresholds, create a plot that displays the Height
signal for the threshold channel and observe the effect on data as the threshold
is raised or lowered.

 Table 3-2.
 Suggested Starting FSC-H Threshold Settings for Various Cell Types

Cell Type	Start with FSC-H Threshold of:	Notes
Cell Lines: Large (>20 µm)	500,000 to 1,000,000	Displaying FSC and SSC signals on Log scale is helpful
Cell Lines: Small (<20 µm)	200,000 to 500,000	Displaying FSC and SSC signals on Log scale is helpful
Unfixed, freshly isolated cells: white blood cells, spleen, thymus	200,000 to 500,000	Linear or log scale display of FSC, SSC
Fixed cell suspensions: white blood cells, spleen, thymus	200,000 to 500,000	Linear or log scale display of FSC, SSC
Platelets	10,000 to 30,000	Linear or log scale display of FSC, SSC
Bacteria, microparticles	Dual thresholds suggested	See document titled "Threshold and Analysis of Small Particles" on the BD Accuri website

**CAUTION:** Take care when setting thresholds before or during data collection. Any event not meeting the threshold criteria will not be acquired or saved. When changes are made to the threshold values after data collection, BD Accuri C6 Software displays a warning message if the new threshold value will result in permanent data loss.



Figure 3-3. Threshold Settings Warning Message

To set the threshold:

- Do one of the following:
  - a. Select Instrument > Set threshold.
  - b. Click on the Set Threshold button in the Instrument Control Panel.

• Select the primary threshold parameter from the *Primary Threshold* drop-down list in the Threshold Settings dialog box.

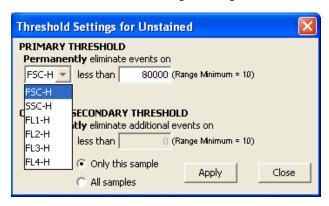


Figure 3-4. Primary Threshold Drop-Down List

• Type 80000 in the less than edit box to set the threshold minimum to channel 80,000.

**NOTE:** A lower or higher FSC-H threshold may be required when working with small cells (such as platelets or bacteria) or large cells (such as cell lines), respectively. Refer to <u>Technical Bulletin: Threshold and Analysis of Small Particles on the BD Accuri C6 Flow Cytometer</u>

(http://www.accuricytometers.com/resources/application-notes/)or Table 3-2.

To apply a secondary threshold:

- Select the threshold parameter from the Secondary Threshold drop-down list.
- Type a value in the less than edit box to set the threshold minimum.
- Do one of the following:
  - a. Select the *Apply to All samples* radio button to apply settings to all samples, including all previously collected data in other data wells.
  - b. Select the *Apply to Only this sample* radio button to apply settings to the current sample only.



Figure 3-5. Threshold Settings Dialog Box

Click on the *Apply* button to apply the threshold settings.

Click on the *Close* button to close the dialog box.

#### 3.2.3 Setting a Run Limit

The run limit defines when data acquisition will stop. The following parameters can be used individually or in combination to set a run limit.

- Time
- Volume
- Number of events (in a specified gate)

If multiple run limits are set, data collection stops on whichever limit is reached first.

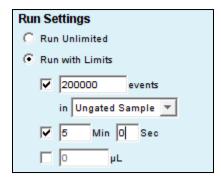


Figure 3-6. Run Limits Controls

To collect samples without setting a run limit:

• Enable the Run Unlimited check box. This requires a manual stop.

To stop the run after a specified number of events have been sampled:

- Enable the Run with Limits check box.
- Enable the check box next to the events field.
- In the associated text box, type the number of events at which to stop the run.
- Do one of the following in the drop-down list below the text box:
  - a. Select Ungated Sample.
  - b. Select a gating strategy (if one exists) to stop the run when the assigned number of events has been collected in the gated region.

To stop the run after a time has expired:

- Enable the Run with Limits check box.
- Enable the check box next to the min and sec fields.
- Type the number of minutes (min) and seconds (sec) at which to stop the run.

To stop the run after a specified volume has been sampled:

- Enable the Run with Limits box
- Enable the check box next to the μL field.
- Type the volume in microliters (µL) at which to stop the run.

#### 3.2.4 Naming the Sample

Samples can be named at any time. If no text is entered in the naming field, BD Accuri C6 Software names the sample according to the well location (for example, A01).

To name the sample:

Type the sample name into the text box above the 96-well grid.



Figure 3-7. Sample Name Field

# 3.2.5 Running the Sample

- Gently resuspend the cells in the sample tube and place the tube on the SIP.
- Select an empty sample well in the BD Accuri C6 Software Collect tab.
- Click on the RUN button to start the sample collection.
  - BD Accuri C6 Software begins fluidics initialization. During this time the Traffic Light turns yellow and the software displays the message *Preparing to analyze sample*. Once initialization is complete, the Traffic Light turns green and the software displays the message *Events are being recorded*. The current well flashes blue during data collection. After the run limit is reached, the well stops flashing and remains blue, indicating that the well contains data.
- Additional data can be collected in a well that contains data by clicking on the ADD TO button. Note that run limits may need to be adjusted when adding data to a well.

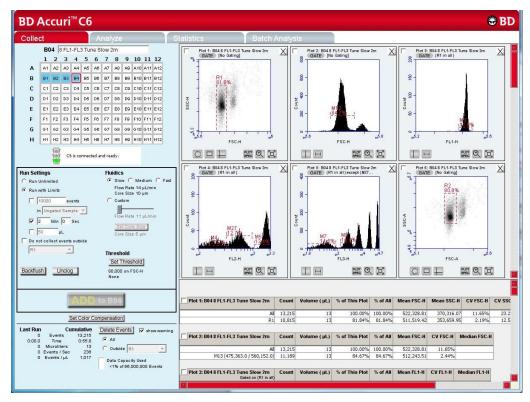


Figure 3-8. BD Accuri C6 Software Workspace after Collecting a Sample

# 3.2.6 Adding More Sample Data to a BD Accuri C6 Software File

Additional sample data can be added to a BD Accuri C6 Software file that already contains data, either by moving to an empty data well before acquisition or by adding to a well which already contains data.

To add data to a BD Accuri C6 Software file:

Gently resuspend the sample and place the tube on the SIP.

**NOTE:** Generally, there is no need to perform a backflush between samples.

Click on a data well in the 96-well sample grid. If an empty well is selected, any
plots and gates created previously are still displayed, but they do not contain any
data, as shown in the following figure.

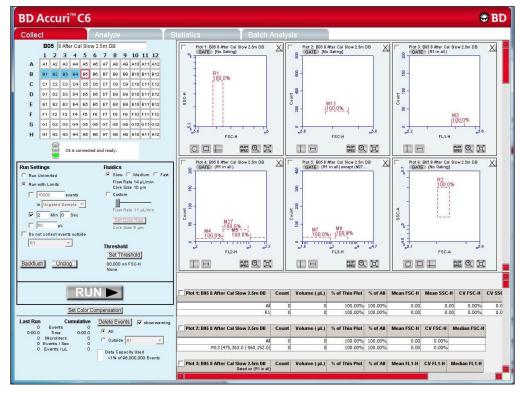


Figure 3-9. BD Accuri C6 Software Workspace with Empty Sample Well and Empty Plots

Click on the RUN (or ADD TO) button to start sample collection. BD Accuri C6
Software displays and updates data in real-time during the collection. The
cytometer stops sampling from the tube when the run limit is reached.

**CAUTION:** If the *ADD TO* button is selected, BD Accuri C6 Software will collect data into a well that already contains data.

### 3.2.7 Pausing Data Collection

Sample acquisition can be interrupted any time during a run. To stop a run:

Click on the PAUSE button.

#### To restart the run:

 Click on the ADD TO button. BD Accuri C6 Software resumes data collection in the current well.

### 3.3 Ending a Data Collection Session

When sample collection is completed, clean the SIP and fluidics lines by following the procedure below:

- Place a tube with 2 mL of filtered, deionized water on the SIP and advance to any empty data well.
- Set the time limit for two minutes and the fluidics speed to Fast.
- Click on the RUN button.
- Place a tube with 2 mL of decontamination solution (PN 653154) (dilute per label instructions) on the SIP.
- Select an empty data well.
- Set the time limit for two minutes and the fluidics speed to fast.
- · Click on the RUN button.
- Once the run is finished, remove the tube of decontamination solution from the SIP.
- Place a tube with 2 mL of filtered, deionized water on the SIP and advance to any empty data well.
- Set the time limit for two minutes and the fluidics speed to Fast.
- Click on the RUN button.
- When the run is finished, leave the tube on the SIP.

### 3.4 Creating Plots

Three plot types are available for viewing data: histogram, density, and dot plots.

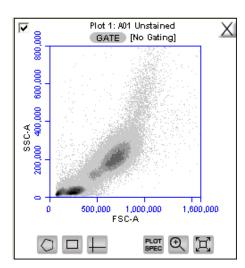


Figure 3-10. View of a Plot

Each plot contains a set of gating and marker tools, and a set of viewing tools:


- Gating and marker tools:
  - Gate button GATE —Opens the Change Gating dialog box for applying gates to a plot.
  - Polygonal Gating Tool For drawing irregularly shaped gates around a population of events.
  - Rectilinear Gating Tool For drawing a rectilinear gate around a population of events.

  - Vertical Marker Tool For gating histograms to the right or left of a vertical marker.
  - Horizontal Marker Tool For gating histograms within a horizontal marker.
- Viewing tools:
  - Plot Spec Tool —Opens the Set Plot Specs dialog box for changing the x- and y-axis parameters, scaling the plot, and setting log or linear view.
  - Zoom Tool —Defines the Zoom range.
  - Expand Tool 
    —Undoes one Zoom level.

#### To create a new plot:

- Click on one of the following icons in an empty plot corral:
  - Density Plot
  - Dot Plot
  - Histogram Plot

BD Accuri C6 Software displays an FSC-A vs. SSC-A plot (or FSC-A, for histogram) by default.

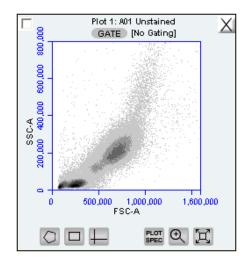

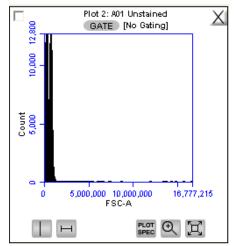


Figure 3-11. New Density and Histogram Plots

Configure the plot specifications as needed.

### 3.5 Changing Plot Specifications

The Plot Spec Tool allows manipulation of the data display in a plot, including axis parameter selection, channel range specification, and selection of linear or logarithmic axis scale. The Plot Spec Tool is available in the *Collect* and *Analyze* tabs.

Set up or modify plot specifications at any time before or after collecting data.

To change the plot specifications:

• Click on the Plot Spec Tool I icon.

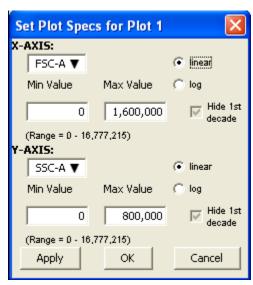


Figure 3-12. Set Plot Specs Dialog Box

- In the Set Plot Specs dialog box, do the following for each axis:
  - Select the desired parameter for display from the drop-down list.
  - Select the *linear* or *log* radio button to specify how data are scaled. (All parameters are collected with linear amplifiers on the cytometer, and therefore all channel values are linear. Selecting the log view of a parameter simply means the linear channel values are graphed on a logarithmic scale.)
  - Type in the minimum and maximum channels to set the channel range to view.
  - Enable or disable the Hide 1st decade check box. It is often beneficial to unhide the first decade of a given parameter before applying fluorescence compensation. See section 3.13.
- Do one of the following:
  - Click on the Apply button to apply the changes without closing the dialog box.
  - Click on the OK button to apply the changes and close the dialog box.
  - Click on the *Cancel* button to close the dialog box without applying the changes.

### 3.6 Using Gates and Markers

A gate is a specified area within a plot that is used to designate a set of events for analysis. BD Accuri C6 Software allows creation of any of the following types of gates:

- Polygonal gate—Gates an irregularly shaped area around a population of events.
- Rectilinear gate—Gates a rectilinear area around a population of events.
- Quadrant gate—Gates the plot in quadrants.
- Vertical marker—Gates a histogram plot to the right or left of a vertical marker.
- Horizontal marker—Gates a histogram plot within the upper and lower boundary of a horizontal marker.

### 3.6.1 Creating a New Gate

To create a gate in a density or dot plot:

- Click on one of the following gating tools:
  - Polygonal Gating Tool typically used for irregularly shaped populations.
  - Rectilinear Gating Tool □—typically used for evenly shaped populations.
  - Quadrant Gating Tool 

    —typically used for analyzing fluorescence plots.
- Use the mouse to draw a region (labeled *P1* for a polygonal gate, *R1* for a rectilinear gate, or *Q1* for a quadrant gate). To draw a polygon, click on the mouse to anchor each vertex and double-click to close the polygon.

**NOTE:** Gate labels can be changed by double-clicking on the label and typing a new gate name in the dialog box.

BD Accuri C6 Software automatically displays the percentage of cells within the region.

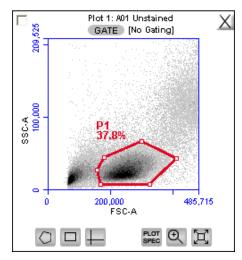


Figure 3-13. Using Polygonal Gating Tool

To create a vertical marker in a histogram plot:

- Click on the Vertical Marker Tool
- Click the cursor at the point along the x-axis on which to place the marker. BD
   Accuri C6 Software automatically displays the percentage of cells to the left (V1-L) and right (V1-R) of the marker.

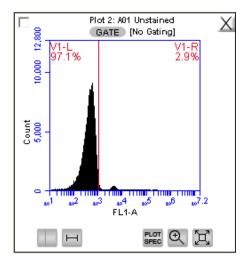


Figure 3-14. Using the Vertical Marker

To create a horizontal marker in a histogram plot:

- Click on the Horizontal Marker Tool .
- Click and drag the cursor horizontally across the area to apply the gate. BD
   Accuri C6 Software automatically displays the percentage of cells within the
   margins of the marker (labeled M1).

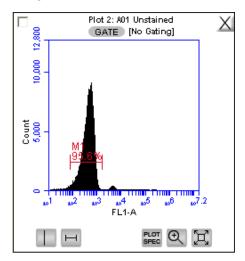


Figure 3-15. Using the Horizontal Marker

# 3.6.2 Applying a Gate to a Plot

- Click on the *GATE* button at the top of the plot. Only polygon (P), rectilinear (R), and marker (M) gating regions automatically appear in the Gating dialog box list of options.
- To include vertical markers or quadrant markers in the list of gates, enable the associated check box(es) in the Change Gating dialog box.



Figure 3-16. Selecting a Gating Option

- Select one of the following gating icons:
  - *Include* icon ——to analyze the events within the region. Multiple gates can be included in a single plot.
  - Exclude icon ——to analyze the events outside of the region.
  - Intersection icon ——to analyze the events within the intersection of two or more regions.
- Click on the Apply button. BD Accuri C6 Software displays the type of gate that is applied next to the GATE button in the plot.

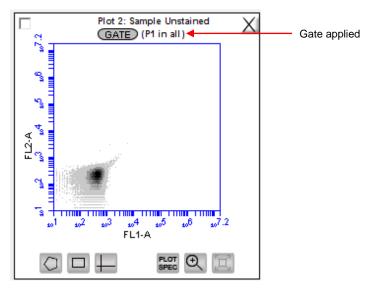


Figure 3-17. Plot Gated to Include P1

# 3.6.3 Creating and Applying Nested Gates

A series of nested gates can be created in which each gate is a subset of the previous one. Complex and informative gating strategies can be devised by the appropriate combination of nested gates using the Include, Exclude and Intersection gating tools.

To create nested gates:

• Draw any region or marker around a population of events (for example, P1).

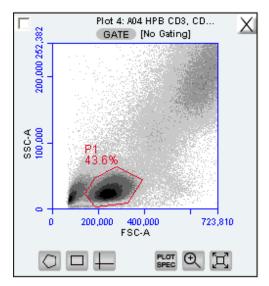


Figure 3-18. First Gate for Creating Nested Gates, Drawn in Plot 4

• Create a new plot and click on the GATE button to apply the P1 (parent) gate

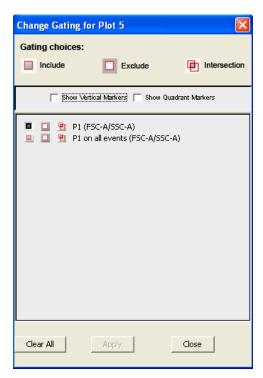


Figure 3-19. Applying the Parent Gate to Plot 5

 Close the dialog box. The plot displays only the populations within the parent gate.

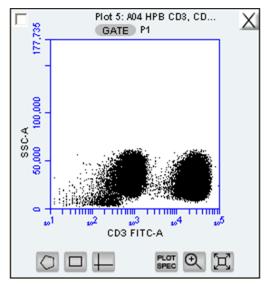


Figure 3-20. Parent Gate Applied to Dot Plot

• In the plot that is gated on the parent gate (plot 5 in this example), draw a second region or marker around a subset of the population displayed in the plot (R1).

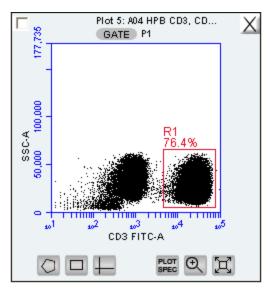


Figure 3-21. Second Gate for Creating Nested Gates

- Create a third plot and click on the GATE button.
- In the Change Gating dialog box, select the option in which the second gate is "in" the parent gate (e.g., R1 in P1; see Figure 3-22). Alternatively, selecting the on all events option (e.g. R1 on all events) "un-nests" the gates.

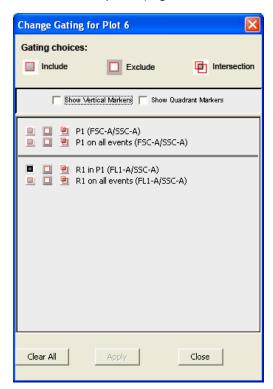


Figure 3-22. Applying the Child Gate

Apply the gate. This is the child gate.

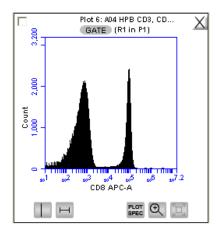


Figure 3-23. Third Plot with Nested Gate Applied (R1 in P1)

Statistics that reflect the nested gates can be viewed in the Statistics Table...



Figure 3-24. Statistics of Plot with Nested Gate

· Close the dialog box.

For instructions on renaming plots and regions, and event coloring see APPENDIX F, "Enhanced Analysis Features."

# 3.7 Moving and Resizing Regions

Region Labels can be moved by clicking and dragging the label. To move or resize a region:

· Click on the border of the region.

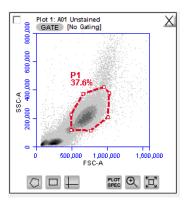


Figure 3-25. Region P1 in the plot has been selected, as indicated by bolded outline

Drag the region according to the desired position.

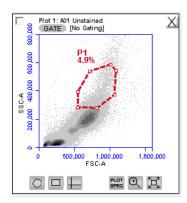
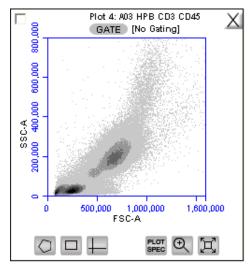


Figure 3-26. Moved region in the Batch Analysis tab

### 3.8 Changing the Number of Events in a Plot

The number of events displayed in all plots across all samples can be changed in order to improve data visualization or to normalize data sets. This option allows visual removal of a number of events from the plot without deleting data.



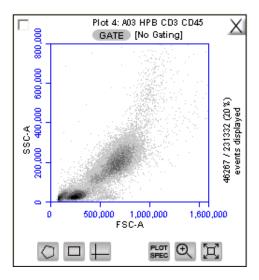


Figure 3-27. Before and After Changing Events Displayed

To change the number of events displayed in a plot:

- Select Display > Events Display Settings.
- Do one of the following in the Events Display Settings dialog box:
  - To view all collected events, select the Show all events radio button.
  - To view the first N events of a sample, select the *Display first* radio button and type a number in the *events collected* field. (Use this option for data normalization.)
  - To view a specified percentage of the whole in a pseudo-random selection, select the *Display* radio button and type a percentage to view (for example, if 20% is selected, every fifth event is displayed).



Figure 3-28. Events Display Settings Dialog Box

- Do one of the following:
  - Click on the Apply button to apply settings without closing the dialog box.
  - Click on the OK button to apply settings and close the dialog box.
  - Click on the Cancel button to close the dialog box without applying settings.

BD Accuri C6 Software displays the plot with a message that some events are not being displayed.

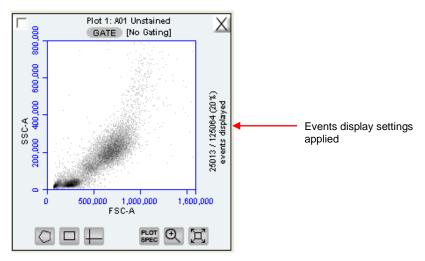


Figure 3-29. Plot with Events Display Settings Applied

# 3.9 Naming Plot Axes

The axes labels can be renamed in any plot from the *Collect* or *Analyze* tab to identify the antibody staining or fluorochrome used in the sample.

To name a plot axis:

• Click on an axis label and select *Rename Parameters* from the drop-down menu. *Rename Parameters* can also be accessed from the *Edit* drop-down menu.

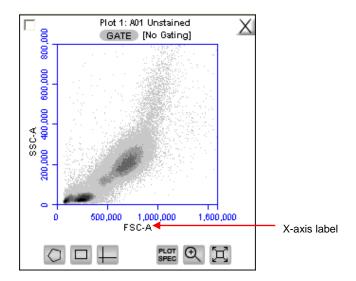


Figure 3-30. X-Axis Label

 In the Rename Parameters dialog box, type the new label in the edit box of the parameter. Any parameter can be renamed from this dialog box.

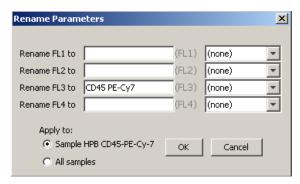


Figure 3-31. Rename Parameters Dialog Box

- Do one of the following:
  - Select the Sample XXX radio button (where XXX refers to the current sample) to apply the label to the current sample only.
  - Select the All Samples radio button to apply the label to all samples.
- Click OK.

To assign the name to the same parameter in another sample:

- Select another sample from the 96-well grid.
- Click on an axis label in a plot and select *Rename Parameters* from the drop-down menu.
- Select the name from the drop-down list associated with the parameter.

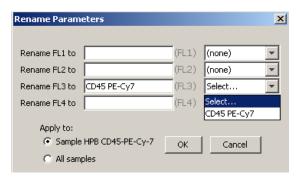


Figure 3-32. Rename Parameters Dialog Box with Axis Label Drop-Down List

#### 3.10 Zooming on a Plot

BD Accuri C6 Software automatically zooms the initial display of any parameter on a logarithmic scale from channel 10 to 16.7 x10<sup>6</sup>. For most analyses, very few events fall into channels 0 to 10, so automatic zooming saves time by reducing the number of zoom steps. However, take care when setting markers (M), regions (R), or polygons (P) that require channels lower than 10 on a zoomed plot. It is recommended to unhide the first decade of data (see section 3.5) to prevent events from being excluded from gated regions, especially when setting fluorescence compensation (see section 3.13).

#### 3.10.1 Basic Zoom

- Click on the Zoom Tool 

  .
- Click and drag the mouse in the plot to draw a zoom area (Figure 3-33).

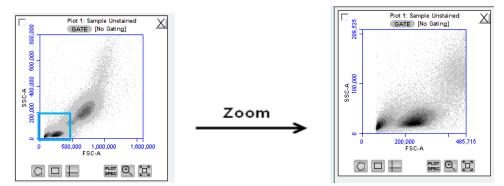


Figure 3-33. Before and After Using Zoom Tool

Repeat steps 1-2 as needed to zoom in.

#### To zoom out:

- Click on the Expand Tool .
- Repeat step 1 as needed.

# 3.10.2 Zooming to a Specified Channel Range

Sometimes it can be helpful to view a plot in a specified channel range.

To view a specified channel range in a plot:

Click on the Plot Spec Tool ....

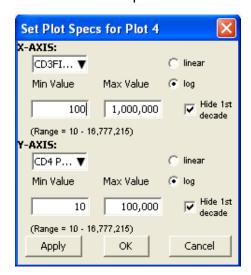


Figure 3-34. Plot Spec Dialog Box: Set Min and Max Channel Values for the X- and Y-Axes

- Specify the x-axis channel range by typing a minimum (*Min Value*) and maximum (*Max Value*) value under *X-Axis* in the Set Plot Specs dialog box.
- Specify the y-axis channel range by typing a minimum and maximum value under *Y-Axis*.
- Click on the Apply button to apply the changes and click on the OK button to close the Plot Spec dialog box.

# 3.11 Saving a BD Accuri C6 Software File

A BD Accuri C6 Software file is a comprehensive (and often large) data file that contains instrument settings, FCS files, and plot layouts.

The file contains the entire BD Accuri C6 Software workspace, including the following elements:

- Sample data
- Plot layouts
- Zoom levels
- Gating
- Color compensation
- Threshold settings
- Collect tab settings
- Changes made in the Analysis or Statistics tabs

By default, BD Accuri C6 Software automatically saves data at the end of each sample run. Data can also be saved manually. To save the entire BD Accuri C6 Software file, save the file manually (see section 3.11.2).

When a BD Accuri C6 Software file is saved, the software displays the file name in the upper left corner of the workspace (Figure 3-35).



Figure 3-35. Title Bar with File Name

#### 3.11.1 Auto-Saving Files

By default, BD Accuri C6 Software automatically saves the event data any time the cytometer reaches a run limit or if a run is paused. <u>Auto-save does not save changes to acquisition settings, plots, or gating strategies that occur after the initial save when naming the file.</u>

**CAUTION:** If changes are made after a run is paused or completed, BD Accuri C6 Software *does not* automatically save the file; save these changes manually (see section 3.11.2).

To enable or disable auto-save:

- Select File > Auto-save Settings.
- Do one of the following in the Auto-save Settings dialog box:
  - Select the Auto-save Enabled radio button to enable auto-save.
  - Select the *Auto-save Disabled* radio button to disable auto-save.



Figure 3-36. Auto-Save Settings Dialog Box

- Click on the OK button to accept the change and close the dialog box.
- If prompted to save the workspace before closing, do one of the following:
  - Click on the Yes button to save the entire workspace.
  - Click on the *No* button to exit the dialog box without saving the workspace.

# 3.11.2 Manually Saving Files

To manually save a BD Accuri C6 Software file:

• Select File > Save.

To manually save a BD Accuri C6 Software file with a new name:

Select File > Save Workspace As.

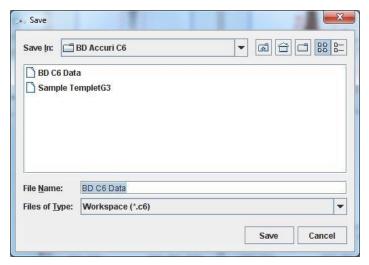


Figure 3-37. Save BD Accuri C6 Software File

- If necessary, navigate to the location to save the file.
- In the Save dialog box, enter the file name and click on the *Save* button. The file is saved with the extension .c6.

# 3.12 Creating a BD Accuri C6 Software Template

A BD Accuri C6 Software template contains a predefined BD Accuri C6 Software workspace for quick and easy setup and analysis. All markers, regions, gates, parameter names, and sample names are saved without any data points. Several templates are provided on the BD Accuri C6 Software CD, flash drive or the BD Accuri website at <a href="www.accuricytometers.com/resources/templates/">www.accuricytometers.com/resources/templates/</a>) or custom templates can be created.

To create a template:

- Define plot, gating, and acquisition settings in a blank workspace, or use the current .c6 file.
- Select File > Save Template as.

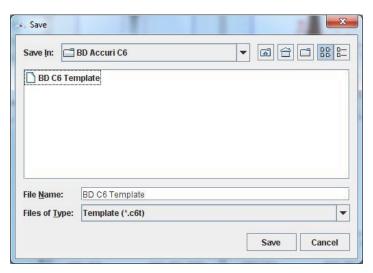


Figure 3-38. Save BD Accuri C6 Software Template

- If necessary, navigate to the location to save the template file.
- In the Save dialog box, enter the file name and click on the Save button. BD Accuri C6 Software saves the file with the extension .c6t.

**NOTE:** Acquisition settings are saved based on the currently selected sample.

### 3.13 Understanding Fluorescence Spillover

Fluorochromes typically emit light over a broad range of wavelengths, resulting in the fluorescence signal appearing not only in the expected, primary detector for that fluorochrome but in other detectors as well. This phenomenon is often called fluorescence "spillover," and can be a source of confusion when interpreting multicolor flow cytometric data.

# 3.13.1 Recognizing Fluorescence Spillover

When performing a multi-color experiment, prepare a set of control samples stained with individual fluorochromes used in the experiment. These single-stained controls will allow determination of the extent of fluorescence spillover from each fluorochrome. The example shown in the figure below shows data collected for a PE-Cy7 single-stained control (red line in Overlay plots) compared to that of unstained cells (black line in Overlays). Most of the fluorescence signal from PE-Cy7 positive cells is detected in the FL3 (670 LP), as expected. However, there is also PE-Cy7 signal detected in FL1 (530 or 533 BP) and FL2 (585 BP), so that plots of data for those detectors appear to have positively fluorescent cells. No signal from PE-Cy7 appears in detector FL4.

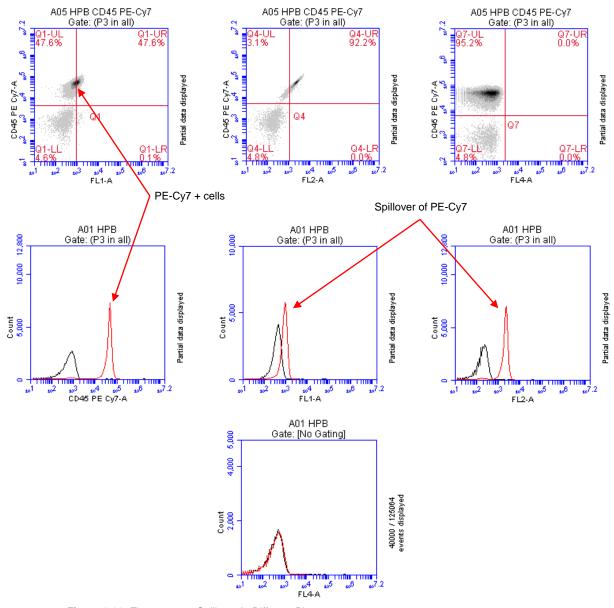


Figure 3-39. Fluorescence Spillover in Different Plots

# 3.13.2 Correcting Fluorescence Spillover

Fluorescence spillover can be removed from plots by applying a mathematical algorithm to collected data. This process is often called *color compensation* (or *fluorescence subtraction*). Because data collection on the cytometer is digital, color compensation can be applied or removed before, during, or after data collection. The color compensation algorithm subtracts a user-defined percentage of fluorescence signal from every event, thereby redistributing data to lower channels on the fluorescence scale and removing the apparent fluorescence spillover. When color compensation has been properly applied to a data set, the median fluorescence channel value in non-primary detectors for any given single-stained control sample should be the same as that of an unstained control sample.

The following figure shows the data after proper color compensation has been applied to the PE-CY7 example in section 3.13.1:

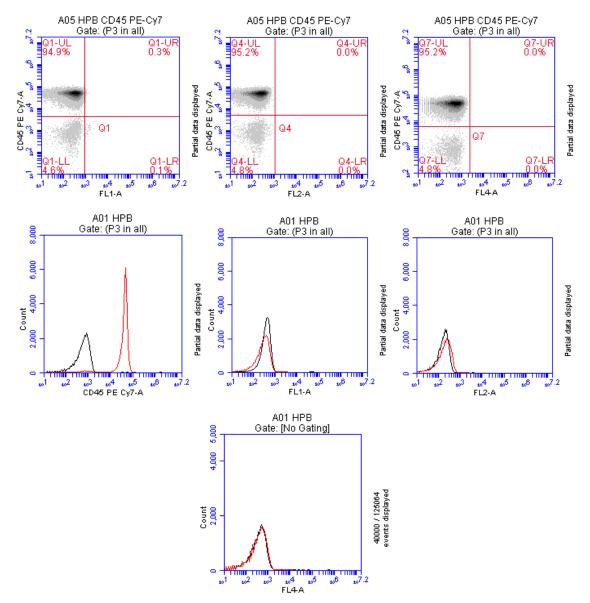


Figure 3-40. Corrected Fluorescence Spillover

The PE-CY7 fluorescence is now confined to the FL3 detector, and no longer spills into FL2 or FL1.

To correct fluorescence spillover:

Click on the Quadrant Tool 

of a plot and click inside the plot.

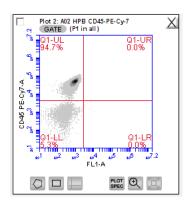


Figure 3-41. Placing a Quadrant Tool

 Adjust the quadrant marker position so that all positive populations are cleanly contained in individual quadrants.

BD Accuri C6 Software displays the median fluorescence channel value for the events in each quadrant in the Statistics Table (shown below). (Be sure that the Show Median Statistics option under the Display dropdown has been selected.)

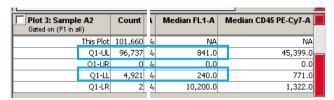


Figure 3-42. Statistics Table Displaying Median Values

- Compare the median values of the affected channel. If the median value of the UL or LR quadrant is not equal to the median value of the negative population (LL), fluorescence compensation should be applied.
- Click on the Set Color Compensation button in the Collect or Analyze tab to open the Compensation Settings dialog box. The dialog box contains four rows of FL buttons, one row for each fluorescence channel.

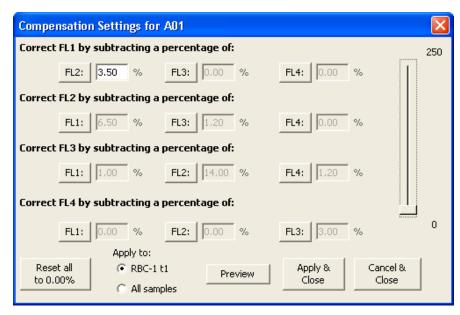


Figure 3-43. Compensation Settings Dialog Box

- In the row associated with the channel to correct, click on the *FL* button of the fluorescence channel that is spilling over.
- If needed, click on the Reset all to 0.00% button to clear all compensation values.
- · Do one of the following:
  - In the text box next to the *FL* button, type an arbitrary percentage of the signal to subtract.
- Use the C Comp Calculator Excel spreadsheet provided on the BD Accuri C6 Software CD (or flash drive) or at <a href="https://www.accuricytometers.com/resources/manuals/">www.accuricytometers.com/resources/manuals/</a>) to calculate the subtraction values.

See the appropriate tab of the C Comp Calculator for your instrument serial number for suggested fluorochrome-specific spillover values (Table 3-3 is an example).

Table 3-3. Example of Fluorochrome Spillover per Channel,

		Spillover Fluorochrome					
		FITC	PE	PerCP	PerCPCy5.5	PE-Cy7	APC
ect	<b>FL1</b> (530BP)	N/A	3.2	0.00	0.0	0.5	N/A
to Correct	<b>FL2</b> (585 BP)	7.5	N/A	0.00	0.00	1.5	0.0
nel t	<b>FL3</b> (670 LP)	1.0	19.5	N/A	N/A	N/A	0.8
Channel to	<b>FL4</b> (675 BP)	0.0	0.0	3.00	12.00	0.00	N/A

- Click on the Preview button to update the plots and Statistics Table.
- In the Statistics Table, observe the values in the Median column of the FL channel of interest.

 Repeat steps 5-9 until the median value for the UL or LR quadrant is equal (or nearly equal) to the median value of the negative population (LL). This value is called the *compensation value*. The figure below shows the median values highlighted in blue.




Figure 3-44. Results of Subtracting Spillover

- To apply the fluorescence subtraction to all samples, select the Apply to All samples radio button in the Compensation Settings dialog box.
- Click on the Apply & Close button to apply the color compensation settings.

### 3.13.3 Troubleshooting Color Compensation

Occasionally, the plot might appear to have a smaller percentage of events in a quadrant than BD Accuri C6 Software reports. For example, Figure 3-45 appears to have less than 26.1% of the population in the lower right quadrant, even though the statistics report 26.1%. This occurs when a decade containing data is hidden.

In the following figure, a number of events have been driven into channel 1 because of overcompensation. These events are not displayed in the plot if the first decade is hidden, but they are included in statistics calculations.

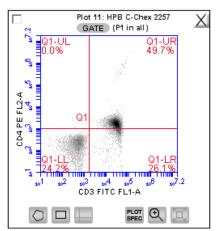

Plot 11: HPB C-Chex 2257 Gated on (P1 in all)	Count	Volume ( μL)	% of This Plot
This Plot	21,479	73	100.00%
Q1-UL	7	73	0.00%
Q1-UR	10,669	73	49.70%
Q1-LL	5,198	73	24.20%
O1-LR	5,605	73	26.10%

Figure 3-45. Plot Displaying Overcompensation


To fix overcompensation:

- Click on the Plot Spec Tool ....
- In the Set Plot Specs dialog box, disable the *Hide 1st decade* check boxes for both X- and Y-axes.



Figure 3-46. Set Plot Specs Dialog Box with Hide 1st decade Disabled

- Click on the OK button to apply the settings and close the dialog box.
- Open the Compensation Settings dialog box and reset the appropriate compensation values to zero. In this example, that would be "Correct FL2 by subtracting a percentage of FL1".
- Perform the fluorescence compensation procedure described in section 3.13.2.
   The number of events displayed in Q1-LR agrees with the percent value of 18.4%, as shown on the plots and Statistics Table in Figure 3-47.

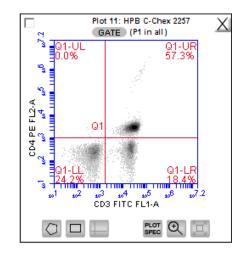

Plot 11: HPB C-Chex 2257 Gated on (P1 in all)	Count	Volume ( µL)	% of This Plot
This Plot	21,479	73	100.00%
Q1-UL	8	73	0.00%
Q1-UR	12,315	73	57.40%
Q1-LL	5,197	73	24.20%
Q1-LR	3,959	73	18.40%

Figure 3-47. Properly Compensated Data after First Decade Is Rehidden

# 3.14 Changing Parameters

By default, BD Accuri C6 Software displays the area parameter (signified with a suffix of -A), but parameters can also be displayed in height, width (of primary threshold channel, only), or time.

To change a parameter:

Click on the x- or y-axis label and select an option from the drop-down menu.

**NOTE:** The time parameter starts counting when the *Run* button is clicked and continues counting for 19 days (16 million tenths of a second), even if data is added to the sample at a later time. The time parameter cannot be reset to zero, even by deleting data.

### 3.15 Copying and Pasting Plots

To copy and paste plots from the *Collect* or *Analyze* tab to a Microsoft<sup>®</sup> Office compatible application:

Click anywhere on a plot and drag it to an open Microsoft application.

**NOTE: Ctrl+C** and **Ctrl+V** cannot be used to copy and paste plots from BD Accuri C6 Software into other applications.

If the Enhanced Analysis activation key is installed, one of two file formats can be selected for plots during drag and drop actions.

To select the file format to use:

format.

- Select File > Set Plot Drag and Drop Format.
- In the dialog box, choose one of the following formats:
   .png when lower resolution is sufficient. Drag and drop to Excel requires.png
  - .eps when higher resolution images are required for publication or posters.



Figure 3-48. Set Plot Drag and Drop Format Dialog Box

**NOTE:** The Drag and Drop Format selection is BD Accuri C6 Software file-specific, and will always revert to the default of .png when a new BD Accuri C6 Software file is created or if the current BD Accuri C6 Software file is not saved with the .eps option selected.

Click on the *OK* button to save the settings and return to the BD Accuri C6 Software Workspace.

### 3.16 Printing Data

To print selected plots and associated statistics from the *Collect* or *Analyze* tab:

Enable the check box in the upper left corner of one or more plots.

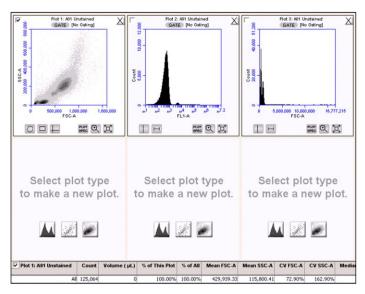


Figure 3-49. Plots Selected for Printing

The associated statistics are automatically selected for printing. This box can be deselected if desired.

Select File > Print Selected Items.

**NOTE:** Printing directly from BD Accuri C6 Software will result in low resolution images even if the .eps option is selected in the Set Plot Drag and Drop Format dialog box.

# 3.17 Exporting and Importing Files

Data can be exported from sample wells from the *Collect* or *Analyze* tab as individual FCS 3.0 files at any time.

**NOTE:** See Appendix F.4 for information about saving plots as publication-quality images.

To export data:

- Do one of the following:
  - Select File > Export FCS File to export and save the currently selected data well as an FCS 3.0 file.
  - Select File > Export ALL Samples as FCS to export and save all data wells as individual FCS 3.0 files.
  - Select File > Export ALL Samples to Third Party to export and save all data wells as individual FCS 3.0 files that enable auto scaling in third party applications (such as FlowJo).
  - Select File > Export Plot Data as CSV to save an individual file in .csv format. (See APPENDIX H for an explanation of .csv format.)

If prompted to confirm the export, click on the OK button.

To import an FCS data file into BD Accuri C6 Software:

- Select an empty data well in a BD Accuri C6 Software file or template. If the currently selected data well contains data, the software will start FCS file import in the first empty data well.
- Select File > Import FCS File.
- Navigate to the location of the file.

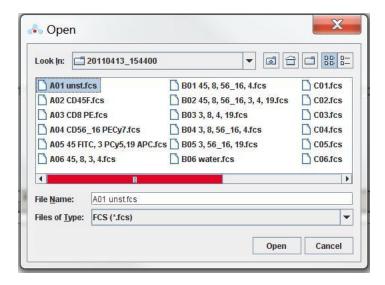


Figure 3-50. Open an FCS file

- Select the file and click on the Open button.
- Multiple FCS files from within a folder can be selected and imported as a group using shift + click. Files begin importing into the selected data well and proceed in succession, horizontally, from left to right, wrapping to the following row once the current row is filled.

**NOTE:** Only FCS files created by BD Accuri C6 Software can be imported into a BD Accuri C6 Software file or template.

### 4 ANALYZING SAMPLE DATA

The *Analyze* tab allows data from multiple samples to be viewed simultaneously using the same plots and gating.

Use the tab to do the following:

- View several plots and samples in any combination for easy analysis.
- Compare specific samples from the 96-well grid.
- Create new plots, hide plots, or copy and reuse plots from the *Collect* tab.
- View different samples with the same plots.
- Create color overlay histograms.
- Print multiple plots.
- Adjust peak position.

# 4.1 Viewing the *Analyze* Tab

The *Analyze* tab is organized into two major sections:

- Setup panel—Panel on the left side of the window that contains controls for selecting samples and plots.
- Data display—Large area on the right side of the window that shows sample data in plots and in a Statistics Table.

When the *Analyze* tab is opened for the first time, the workspace is empty. Plots can be copied from the *Collect* tab or created from scratch. Gating strategies that were set up in the *Collect* tab can be applied in *Analyze*, or new gates can be drawn and new gating strategies set up in *Analyze*.

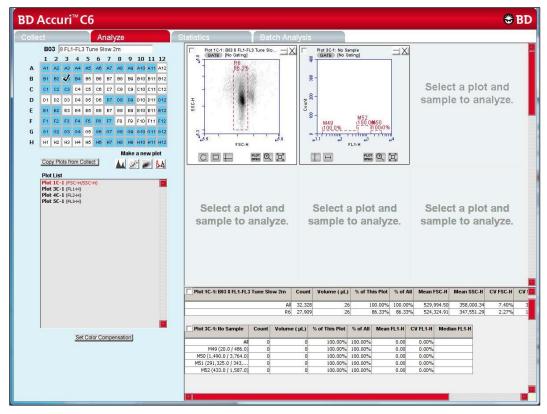


Figure 4-1. Analyze Tab Workspace

The following table describes each of the controls and indicators in the *Analyze* tab:

Table 4-1. Analyze Tab Controls

Control	Description	
Sample Naming Field	Text box for naming the current sample.	
Sample Grid	Matrix laid out in the configuration of a 96-well plate to organize sample data. Each sample has its own well in the Sample Grid.	
	The wells are color-coded:	
	White—Does not contain data.	
	Blue—Contains data.	
	Black check mark—Currently selected for viewing data.	
Copy Plots from Collect	Copies specified plots from the <i>Collect</i> tab. See section 4.2.1 for details.	
Plot Controls	Set of buttons for creating new plots or overlaying histograms. See section 4.2.3 for details.	
	All plots created in the <i>Analyze</i> and <i>Collect</i> tabs, including overlays, can be dragged and dropped into most Microsoft Office compatible programs.	
Plot List	Lists the plots that are available in the <i>Analyze</i> tab.  Available plots include plots copied from the <i>Collect</i> tab or created in the <i>Analyze</i> tab.	

Control	Description
Set Color Compensation	Opens the Color Compensation Matrix for correcting fluorescence spillover. See section 3.13 for details.
Plot Corrals	Area displaying two rows of plot corrals. Scroll up or down to view more plots. For information on creating plots, see section 4.2.2.
Statistics Table	Table below the plots that displays statistical information on individual plots.
	Statistics Tables can be copied into most Microsoft Office compatible programs.

### 4.2 Setting up Plots

In the *Analyze* workspace, plots can be created or copied from the *Collect* tab. Plots that are copied from the *Collect* tab are appended with a "C" (for example, *Plot 1C*).

NOTE: See APPENDIX F.4 for details on saving plots as publication quality images.

### 4.2.1 Copying Plots from the *Collect* Tab

- In the Analyze tab, click on the Copy Plots from Collect button.
- In the Copy Plots from Collect dialog box, do one of the following:
  - Select plots to copy by checking the box(es).
  - Enable the All Plots check box to copy all plots from the Collect tab.

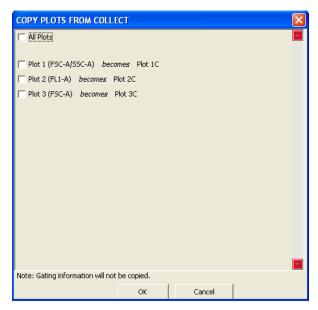


Figure 4-2. Selecting Plots to Copy from the Collect Tab

• Click on the *OK* button to close the dialog box. BD Accuri C6 Software adds the selected plots to the Plot List in the *Analyze* tab (Figure 4-3).

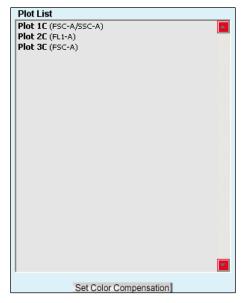


Figure 4-3. Plot List Containing Copied Plots

# 4.2.2 Creating Plots

- Click on an empty plot corral.
- Click on one of the following icons under the Sample Grid:
  - Histogram ▲
  - Dot <u>#</u>
  - Density
  - Overlay Histogram Ы (see section 4.2.3 for details)
- · Click on the sample well to view data.

# 4.2.3 Creating an Overlay Histogram

Creating an overlay histogram allows comparison of multiple distributions, up to 96 different samples at the same time. To create an overlay histogram:

Click on an empty plot corral.

 Click on the Overlay Histogram Tool to open a blank single-parameter FSC-A plot (Figure 4-4).

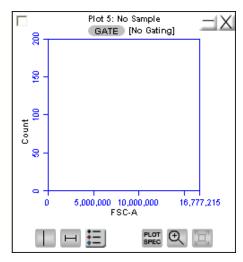


Figure 4-4. Blank Overlay Histogram Plot

- Click on the x-axis label (FSC-A) and select a different parameter in the drop-down list, if desired.
- Click on the GATE button and apply a gate, if desired (see section 3.6 for details).
- Select the samples to be overlaid from the 96-well grid.

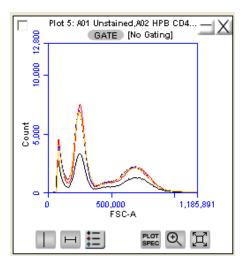


Figure 4-5. Overlay Histogram Plot with Data

Click on the Overlay Histogram Legend Tool 

 is to view a legend for the overlay histogram.

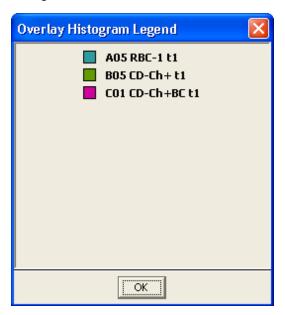


Figure 4-6. Overlay Histogram Legend

To change the color of one or more overlay plots, click on the square in the Overlay Histogram Legend and select the desired color from the pop-up color palette.

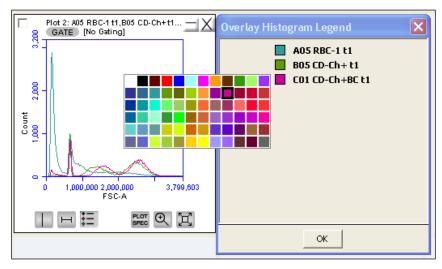


Figure 4-7. Overlay Histogram Legend with Color Palette

# 4.3 Viewing Plots

To view a plot in the Plot List:

Click on an empty plot corral in the Analyze tab.

• Select a plot in the Plot List (Figure 4-8). BD Accuri C6 Software displays the plot without any sample data.

**NOTE:** Any gating tools (markers, regions, etc.) carried over from the *Collect* tab are renamed (for example, P1 in *Collect* is P2 in *Analyze*). When these "new" markers or regions are adjusted in the *Analyze* tab the original marker or region position in *Collect* does not change.

- Click on a sample well to view data.
- Apply a gating strategy, if desired (see section 3.6.2 for details).

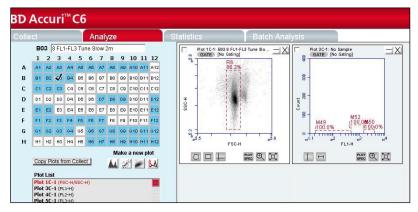


Figure 4-8. Analyze Tab: Plots with Gating Applied

To view data from another sample, open one or more plots from the Plot List (it is recommended to do this in another row of plot corrals) and choose the sample to be displayed in each plot. Gating strategies applied above are automatically applied to the corresponding plots in the new row.

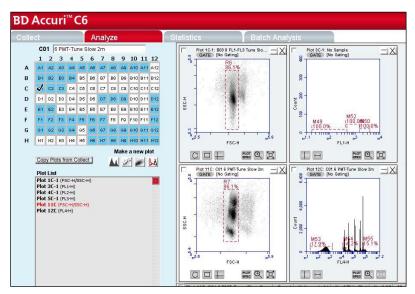


Figure 4-9. Analyze Tab: Two Samples Displayed with the Same Plots

Compare data and statistics between samples.

# 5 CREATING A STATISTICS TABLE

The *Statistics* tab provides a way of tabulating data from multiple samples in one master table. It also allows users to do the following:

- View statistics for some or all samples.
- Display statistics of collected or imported samples.
- List all plots created on the Collect and Analyze tabs.
- Display all plot names, gates, and associated statistics.
- Copy and paste data into a spreadsheet.

### 5.1 Viewing the Statistics Tab

The Statistics tab is organized into two major sections:

- Setup controls for the Master Statistics Table
- Master Statistics Table

When the *Statistics* tab is first opened, the workspace is empty. To create a Master Statistics Table select data from the setup controls.

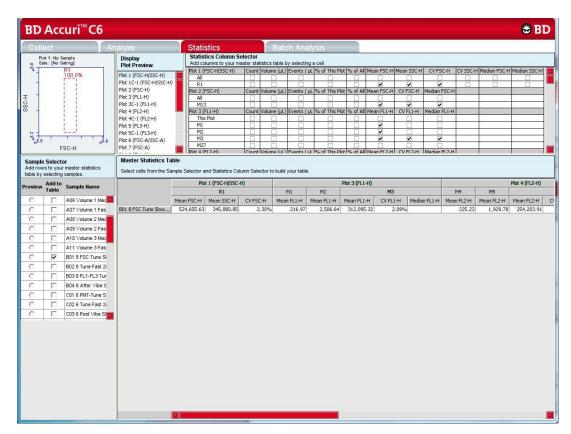


Figure 5-1. Statistics Tab Workspace

The following table describes each of the controls and indicators in the *Statistics* tab:

Table 5-1. Statistics Tab Controls

Control	Description
Plot	Previews the selected plot. See section 5.3 for details.
Display Plot Preview	List of plots imported from the <i>Collect</i> and <i>Analyze</i> tabs that are available to preview.
Statistics Column Selector	Selects the data to view in the Master Statistics Table for each sample. See section 5.2 for details.
Sample Selector	Selects the samples to view in the Master Statistics Table. See section 5.2 for details.
Master Statistics Table	Configurable table that displays data of selected samples. See section 5.2 for details.

### 5.2 Creating the Master Statistics Table

The Master Statistics Table enables creation of a customized data table for multiple samples within a given BD Accuri C6 Software file.

To create the Master Statistics Table:

 In the Statistics Column Selector, enable the check boxes under the data items to view per plot. BD Accuri C6 Software automatically adds columns to the Master Statistics Table.

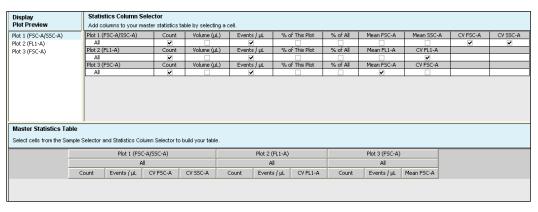


Figure 5-2. Creating Master Statistics Table: Adding Plots

In the Sample Selector list, enable the check box of each sample to be viewed.
 BD Accuri C6 Software automatically adds rows of samples to the Master Statistics Table and displays the sample data.

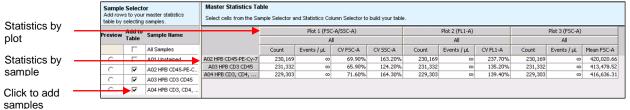


Figure 5-3. Creating Master Statistics Table: Adding Samples

# 5.3 Previewing a Plot in the *Statistics* Tab

To preview a plot in the Statistics tab:

In the Display Plot Preview list, click on the plot to preview.

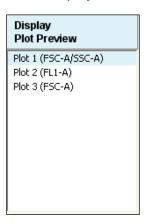


Figure 5-4. Plot Preview

• In the Sample Selector list, select the radio button of a sample. BD Accuri C6 Software displays the sample data in the plot.

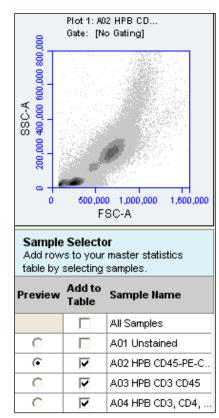


Figure 5-5. Sample Selector List and Plot

The plot preview is available for viewing only. The zoom level and other plot settings must be modified in either the *Analyze* or *Collect* tab.

# 5.4 Copying Data into Other Applications

Data can be copied and pasted from the Master Statistics Table into most Microsoft Office compatible applications.

To copy data:

- Use the mouse to highlight the fields. The column and row headers corresponding to the selected data fields are automatically copied.
- Press Ctrl+C to copy the data.
- In the Microsoft application, press Ctrl+V to paste the data.

**NOTE:** See APPENDIX F.5 for instructions on Batch Analysis.

# 6 MAINTAINING THE BD ACCURI C6 FLOW CYTOMETER

To maintain optimal performance, follow the routines in this chapter on a regular basis. See the *BD Accuri C6 Flow Cytometer Instrument Manual* for mechanical maintenance procedures (such as replacing tubing).

## 6.1 Cleaning the SIP

Run the Backflush Cycle to clean the SIP and remove clogs at the base of the SIP.

To perform a backflush:

- Place a blotter or empty sample tube under the SIP to catch dripping fluid.
- · Do one of the following:
  - Click on the Backflush button in the Collect tab.
  - Select Instrument > Run Backflush Cycle.
- When the backflush is completed, remove the tube from the SIP.

## 6.2 Cleaning the Flow Cell

Clean the flow cell as a part of regular maintenance or to correct performance issues of the cytometer.

# 6.2.1 Running the Unclog Cycle

The Unclog Cycle purges the flow cell of debris.

To purge the flow cell:

- Remove the sample tube from the SIP and place a blotter or empty sample tube under the SIP to catch dripping fluid.
- Do one of the following:
  - Click on the *Unclog* button in the *Collect* tab.
  - Select Instrument > Run unclog cycle.

# 6.2.2 Running an Extended Clean of the Flow Cell

During the Extended Flow Cell Cleaning Cycle, the flow cell fills completely with the extended flow cell cleaning solution from a sample tube on the SIP. This cycle automatically shuts down the cytometer with the solution in the flow cell, allowing the flow cell to soak. It is recommended to perform this cleaning cycle monthly, or when a partial blockage of the SIP or flow cell is suspected.

To run the Extended Clean of the Flow Cell Cycle:

 Place a tube with at least 500 μL of Extended Cell Flow Clean cleaning solution (PN 653159—NOT cleaning solution, PN 653157) on the SIP.

**CAUTION:** Never run the Extended Clean of the Flow Cell Cycle without a tube containing at least 500 µL of fluid.

- Select Instrument > Extended clean of flow cell.
- After the cytometer is shut down, leave the cytometer off for at least 30 minutes (up to overnight, for a more thorough cleaning).
- Restart the cytometer. The cytometer performs a longer fluidics startup cycle and BD Accuri C6 Software displays the message *Extra startup time needed due to cleaning or improper shutdown*. This longer cycle purges cleaning solution from the flow cell and takes about 15 minutes to complete.
- Replace the tube containing cleaning solution with one containing 0.2 µm filtered deionized water, and run on Fast fluidics speed for at least 5 minutes to clear residual cleaner solution from the SIP and flow cell.
- Operate the cytometer as usual.

# 6.3 Cleaning the Fluidics Lines

The Cleaning Fluid Cycle pulls cleaning solution from the cleaner tank and runs it through the fluidic lines. After filling the system with cleaning solution, the cleaning fluid cycle purges the cytometer with fresh sheath fluid and performs a backflush. This cycle takes about five minutes.

To run a cleaning fluid cycle:

- Place a tube of cleaning solution (PN 653157) on the SIP.
- Select Instrument > Run cleaning fluid cycle.

# 6.4 Decontaminating the Fluidics System

BD Accuri C6 Software automatically decontaminates the cytometer fluidics system at shut-down. The Decontamination Fluid Cycle can also be run manually at any time. During this process, the cytometer pulls decontamination fluid from the decontamination bottle then pulls sheath fluid from the sheath fluid bottle. Decontamination takes about 13 minutes.

To manually decontaminate the fluidics:

- Place a tube of water on the SIP.
- Select Instrument > Run decontamination fluid cycle.

# 6.5 Using the BD Accuri C6 Flow Cytometer for Precise Volume Measurements

Each BD Accuri C6 flow cytometer is calibrated at the factory for accurate volume measurement before shipping. Follow the guidelines below when accurate volume measurement is required during sample acquisition. For more detailed information on this subject, refer to the Technical Note, "A Guide to Absolute Cell Counting on the BD Accuri C6 Flow Cytometer" available at www.AccuriCytometers.com.

- Ensure that the peristaltic pump tubing and in-line sheath filter have been replaced within 60 days.
- Use 12x75 mm tubes (any type of plastic). Other tubes are not supported.

- Use sample volumes between 300 μL and 2 mL. Never acquire more than 750 μL from a single tube on *Medium* fluidics setting or 1500 μL on *Fast* fluidics setting.
- Adjust sample concentrations to fall approximately within the range of 1x10<sup>3</sup> and 5 x 10<sup>6</sup> cells or particles per mL.
- Acquire data on Medium or Fast fluidics settings only. Custom fluidics settings above 15 μL and 16 μm may be used, but must be validated by an independent control/count bead.
- Only acquire once from any sample tube. Sample height within the tube is critical. Replicate measurements must be obtained from separate sample tubes made by aliquoting sample into the appropriate number of equal volumes.
- Always compare the same stop limit types. Do not compare concentrations collected with volume stop counts to event stop counts.
- Always validate accurate counting by using a reference count bead in the
  experimental buffer, using the same sample volume and tube as in the
  experiment. If bead counts are within 20% of the expected value (based on
  information provided by the bead manufacturer), proceed with sample collection.
  If bead counts are not within 20% of expected values, proceed with fluidics
  calibration as described below.
- Perform the following, in order:
  - 1. Ensure that the fluid levels in the Sheath, Cleaner and Decontamination bottles are sufficient to cover the inlet tubing and that there are no "kinked" fluidic lines.
  - 2. Run a decontamination cycle from the Instrument Menu or by restarting the cytometer.
  - 3. Within 5 minutes of completion of the water run, place a calibration sample on the SIP. Select Instrument > *Calibrate Fluidics*.
    - Calibration should be performed in the same tube as the experimental sample.
    - Calibration should be performed using a sample of the same or similar viscosity as the samples to be analyzed. For example, if lysed human peripheral blood samples are to be acquired, lysed human peripheral blood should be used during calibration.
    - The volume in the control sample tube should be 110 μL more than the volume used with subsequent test samples. For example, if using 1000 μL samples, perform calibration with 1110 μL in the tube. The calibration procedure consumes approximately 220 μL. The values determined by the BD Accuri C6 are based on the average sample height in the tube during the calibration.
    - If sample volumes >50 μL are to be acquired from the sample tube, the calibration volume should take this into account and the average volume in the sample tube during the acquisition should be used. For example, if 100 μL are to be acquired from a 1000 μL sample, the calibration volume would be 950 μL. The average volume [(Starting Volume + Ending Volume)/2] during the acquisition is [(1000+900)/2].
  - 4. The cytometer will perform a calibration cycle, taking approximately 13 minutes. Once completed, the traffic light will revert to green with the *C6 is*

connected and ready status message. Repeat fluidics validation using a reference count bead as described above.

If calibration fails, BD Accuri C6 Software will display a message that indicates that instrument calibration has failed.

Troubleshooting Instrument Calibration:

- Make sure that the calibration tube did not run dry during calibration.
- Repeat calibration preparation routine, including running the decontamination cycle and water.
- Repeat the calibration with a new calibration sample.
- If calibration fails a second time, replace the peristaltic pump tubing in the cytometer and repeat the preparation/calibration routine.
- If calibration fails a third time, contact BD Accuri Technical Support.
- 5. The cytometer will still operate normally after a failed calibration. However, the volume measurements for the desired samples may be incorrect because the cytometer will revert to the factory-set, default fluidics calibration settings. All other aspects of the data will be normal.

# 7 ADJUSTING PEAK POSITION WITH VIRTUALGAIN™

In certain instances, a particular peak should have the same position across different samples or be located at a specific channel number, regardless of the staining. Instruments that have voltage and amp gain controls allow peak position adjustments from sample to sample. BD Accuri C6 Software uses VirtualGain instead of these controls.

VirtualGain is a software module that mimics voltage and amp gain adjustments to reposition data on the axis after the data has been collected. VirtualGain makes gross adjustments (approximate visual shifts of the data) of histogram plots. It is strictly an analysis tool and should not be used while collecting data.

For example, in Figure 7-1 the negative peaks in the control sample and in sample 1 fall in similar channels (mean value = 28.2 and 29.7, respectively). However, the negative population in sample 2 is farther to the right (mean value = 73.4). VirtualGain can be used to align the negative peak of sample 2 with the control sample.

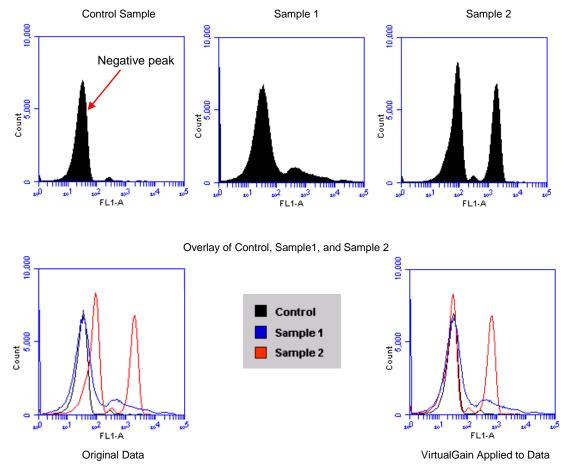


Figure 7-1. Before and After Applying VirtualGain

# 7.1 Applying VirtualGain

Only apply VirtualGain on a histogram plot and only on one parameter at a time. After VirtualGain is applied, data can be viewed in any type of plot and VirtualGain can be toggled on and off.

VirtualGain is only applied to the displayed data and does not alter FCS data. The adjustment is recorded only in the BD Accuri C6 Software file.

### To apply VirtualGain:

- In the Analyze tab, do one of the following:
  - Recreate the histogram to which VirtualGain is to be applied (see section 4.2.3 for details).
  - Copy plots from the Collect tab (see section 4.2.1 for details).

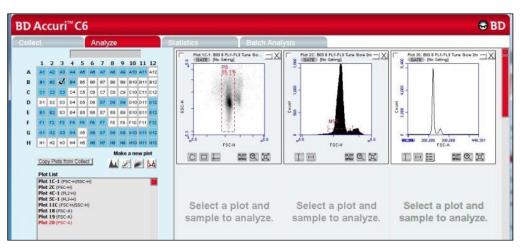


Figure 7-2. Setting Up Histograms for VirtualGain

- Apply the appropriate gating to the plots in the Analyze tab.
- Do one of the following:
  - Select a histogram plot from the sample to which the other samples will be aligned. This sample is the *standard sample*.
  - Select an empty well to align data to a specific channel instead of a collected sample.
- Click on the x-axis label on the standard sample and select *VirtualGain* from the pop-up Parameter List.
- In the VirtualGain dialog box, do one of the following:
  - Move the peak definition marker (vertical line) in the Standard Sample plot to the center of the peak that will be the reference point. Other samples will be aligned to this position.
  - If an empty well was selected in step 3, move the peak definition marker (vertical line) to a channel to assign as a reference point.

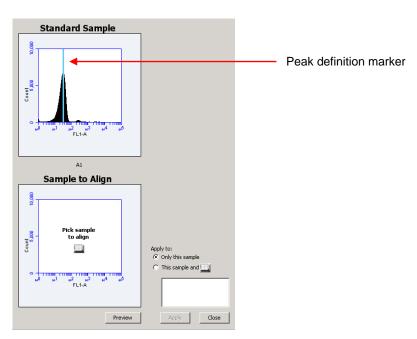


Figure 7-3. Aligning Plots

- If needed, use the Zoom Tools in the *Analyze* tab to change the zoom level in the VirtualGain dialog box.
- Click on the small sample grid icon in the center of the Sample to Align plot.
- Open the sample to be aligned by clicking on the corresponding blue well in the pop-up sample grid (the gray well indicates the standard sample currently selected). Ensure that this plot has been zoomed to the required level before setting VirtualGain.

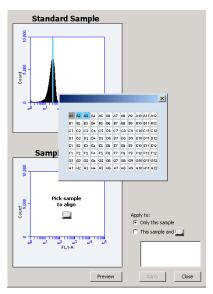


Figure 7-4. Pick the Sample to Align

• Move the peak definition marker in the Sample to Align plot to the center of the peak to align (Figure 7-5).

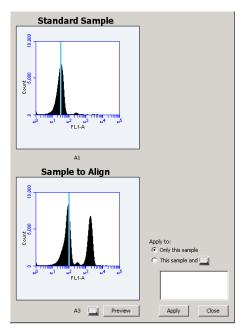


Figure 7-5. Move the Peak Definition Marker

- Click on the *Preview* button to view the aligned sample with VirtualGain applied. BD Accuri C6 Software aligns the peak of interest in both plots.
- Repeat steps 9-10 to make additional adjustments, if needed.
- To align additional samples exactly as the first aligned sample, select the *This* sample and radio button and click on the well(s) in the pop-up sample grid to be included. If the other samples need a different amount of VirtualGain, set VirtualGain separately for each sample.

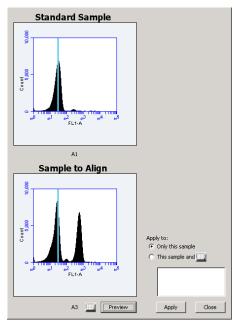


Figure 7-6. Align Additional Samples

 Click on the Apply button to apply VirtualGain to the data. BD Accuri C6 Software displays a black asterisk under the Sample to Align plot to indicate that VirtualGain has been applied to the specified parameter for that sample.

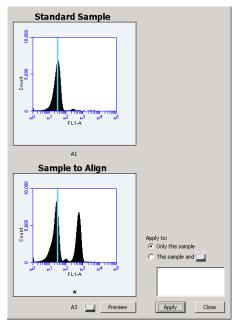


Figure 7-7. Black Asterisk Identifier

• Click on the Close button to close the VirtualGain dialog box.

# 7.2 Viewing VirtualGain

When VirtualGain is applied to a sample, BD Accuri C6 Software displays a black asterisk under the parameter label in the associated plot.

The asterisk is color-coded:

- Black—VirtualGain has been applied.
- Gray—BD Accuri C6 Software is currently displaying the original data.

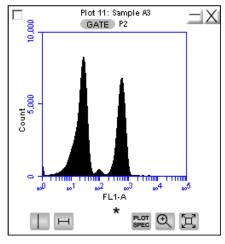


Figure 7-8. Plot with Black Asterisk

Overlays automatically display VirtualGain when it is applied (Figure 7-9). The asterisk is not displayed in overlays when VirtualGain is applied to some or all of the samples in the overlay.

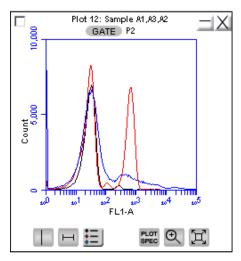


Figure 7-9. VirtualGain Applied in an Overlay Histogram

To toggle between views with VirtualGain applied and not applied:

Click on the asterisk in the plot.

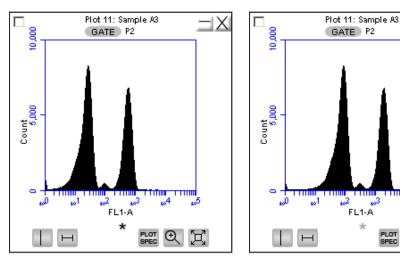


Figure 7-10. Toggle between VirtualGain Applied (Left) and Not Applied (Right)

# 7.3 Removing VirtualGain

To permanently remove VirtualGain from every parameter in every sample in the BD Accuri C6 Software file:

PLOT ① □

Select Display > Remove All VirtualGain.

**CAUTION:** This is a permanent action: Undo cannot be applied.

# APPENDIX A BD ACCURI C6 SOFTWARE MENU QUICK REFERENCE

The following table provides a description of all menu options in BD Accuri C6 Software.

Table A-1. BD Accuri C6 Software Menus

	Accuit Co Software Merius	
Menu Item	Menu pick	Description
File	Open BD Accuri C6 Software File or Template	Opens a previously saved BD Accuri C6 Software file or template. Only one BD Accuri C6 Software workspace can be open at a time.
	New BD Accuri C6 Software File	Opens a new, blank BD Accuri C6 Software workspace. Replaces any previously open workspace.
	Save	Saves the open BD Accuri C6 Software workspace under the current name. If the file has not already been named, BD Accuri C6 Software prompts the user to name the file.
	Save BD Accuri C6 Software File as	Saves the open BD Accuri C6 Software workspace under a new name.
	Save BD Accuri C6 Software Template as	Creates a template from the currently open BD Accuri C6 Software workspace. All markers, regions, gates, parameter names, and sample names are saved without any data points.
	Auto-save Settings	Allows auto-save feature to be enabled or disabled.
	Import FCS File	Imports an FCS file previously exported from another BD Accuri C6 Software file to the currently open workspace. Only FCS files created on a BD Accuri C6 flow cytometer can be imported into BD Accuri C6 Software.
	Export FCS File	Exports and saves the currently selected data well as an FCS 3.0 file to a specified folder. Exported files are compatible with off-line analysis programs such as FCS Express, FlowJo, and WinList.
	Export ALL Samples as FCS	Saves all of the data wells as individual FCS 3.0 files in the folder FCS Exports on the computer desktop. Exported files are compatible with off-line analysis programs such as FCS Express, FlowJo, and WinList.
	Export ALL Samples to Third Party	Exports and saves all data wells as individual FCS 3.0 files that enable autoscaling in third party applications (such as FlowJo).
	Export Plot Data as CSV	Saves an individual file in .csv format for further analysis in spreadsheet programs. All data for every event in the selected plot is exported. See APPENDIX H for an example .csv file.

Menu Item	Menu pick	Description
	Set Plot Drag and Drop Format	Available with Enhanced Analysis Features. Allows plot format to be toggled between .png (low resolution) or .eps (high resolution).
	Print Selected Items	Prints selected plots and associated statistics.
	Quit	Quits BD Accuri C6 Software and closes the application.
Edit	Undo	Undoes the last action that was performed in BD Accuri C6 Software. Not all actions are undoable.
	Redo	Reverses an <i>Undo</i> action.
	Сору	Copies a selected marker or region from a plot, or selected statistics from the tables in BD Accuri C6 Software.
	Paste	Pastes copied markers and regions into new plots.
	Rename Parameters	Allows individual parameters in either the current sample or all samples to be renamed.
Display	Events Display Settings	Opens a dialog to change the number of events displayed in all plots.
	Auto-Select Next Well	Opens a dialog to configure whether BD Accuri C6 Software automatically selects wells vertically, horizontally, or not at all at the completion of each sample collection.
	Remove All VirtualGain	Removes all VirtualGain settings from the entire BD Accuri C6 Software workspace.
	Hide/Show Median Statistics	Hides/shows the median statistics in the Statistics Table.
Instrument	Set threshold	Opens the Threshold dialog box for setting the primary threshold value, and setting an optional secondary threshold.
	Set compensation	Opens the Compensation Settings dialog box for correcting fluorescence spillover.
	Run cleaning fluid cycle	Runs a cleaning fluid cycle.
	Run decontamination fluid cycle	Runs decontamination fluid cycle.
	Run unclog cycle	Runs the unclog cycle to clean the flow cell.
	Run backflush cycle	Runs the backflush cycle to clean the SIP and remove clogs at the base of the SIP.
	Extended clean of flow cell	Cleans the flow cell for an extended time.
	Calibrate fluidics	Initiates fluidics calibration to ensure that the cytometer provides accurate volume measurement.
	Update firmware	Updates the cytometer firmware. Use only when directed to upgrade the cytometer firmware with an official firmware release.
	Remote Control	Optional feature that allows control of the cytometer from a remote location.
About	About Accuri C6 Software	Opens a dialog box that displays the version of BD Accuri C6 Software and BD Accuri Technical Support contact information.

Menu Item	Menu pick	Description
	Technical Support Information	Opens a dialog box that displays information about BD Accuri C6 Software and the BD Accuri C6 flow cytometer. Each time an activation key is used to install a new BD Accuri C6 Software component, the dialog box is updated to reflect the change.
	Users	Allows addition, deletion or modification of user accounts with the optional user tracking feature.

# APPENDIX B ADVANCED FLUIDICS SETTINGS

Advanced users can customize the fluidics rate and core size for collecting samples.

To customize the fluidics rate:

Select the Custom radio button in the Fluidics section of the Collect tab.

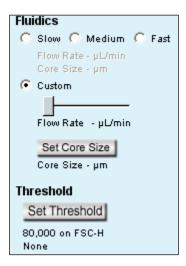


Figure B-1. Custom Radio Button

Move the Custom slider to adjust the flow rate.

To customize the sample core size:

- Click on the Set Core Size button in the Fluidics section of the Collect tab.
- Move the slider to adjust the core size.

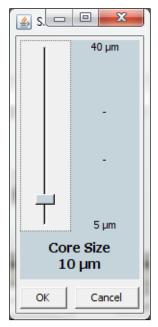


Figure B-2. Set Core Size

**NOTE:** Certain core sizes are not compatible with certain flow rates. BD Accuri C6 Software does not allow these combinations to be set. Use the following table to determine allowable combinations.

 Table B-1. Core Size/Flow Rate Combinations

Core Size	Min Flow Rate	Max Flow Rate
5	10	11
6	10	16
7	10	22
8	10	29
9	10	36
10	10	45
11	10	54
12	10	65
13	10	76
14	12	88
15	14	100
16	15	100
17	17	100
18	19	100
19	22	100
20	24	100
21	26	100
22	29	100
23	32	100
24	35	100
25	38	100
26	41	100
27	44	100
28	47	100
29	50	100
30	54	100
31	58	100
32	61	100
33	65	100
34	69	100
35	74	100
36	78	100
37	82	100
38	87	100
39	91	100
40	96	100

• Click on the OK button to set the core size and close the slider.

# APPENDIX C TRACKING USER ACTIVITY

User tracking allows laboratory administrators to track the activities of BD Accuri C6 flow cytometer operators by assigning a user name and password to each individual. Passwords are created and used in BD Accuri C6 Software and are unrelated to any Windows passwords used on the host computer or network.

User Tracking is an optional upgrade for BD Accuri C6 Software and BD CSampler Software and requires the use of BD Accuri C6 Software version 227 or above.

# **C.1** Installing the User Tracking Module

- Ensure BD Accuri C6 Software version 227 or above is loaded on the computer.
- Copy the User Tracking Activation Key from the installation CD or flash drive to the computer desktop.
- Double-click on the User Tracking Installer icon on the desktop.



Figure C-1. User Tracking Installer Icon

- In the installation wizard, choose the directory in which to install the Activation Key. The location depends on where BD Accuri C6 Software has been installed. In most cases, the location is in
  - C:\Program Files\BD Accuri\BD Accuri C6 Software\ActivationKeys
  - If necessary, use the browse button .... to navigate to the correct folder.

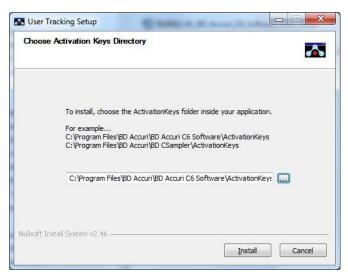


Figure C-2. Installation Dialog Box

 Click on the *Install* button. BD Accuri C6 Software displays a confirmation message after successful installation.

# **C.2** Using the Tracking Feature for the First Time

To use the tracking feature for the first time:

- Open BD Accuri C6 Software.
- When prompted, type admin in the Username text box.

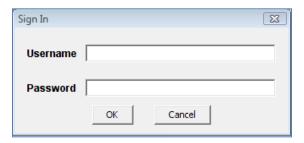


Figure C-3. Username and Password Dialog Box

- Type Admin in the Password text box (case sensitive). The Administrator password can be changed later.
- Click on the OK button
- Use BD Accuri C6 Software as usual.

## C.3 Adding, Deleting, and Modifying User Accounts

The administrator can add new users, delete existing users, and change passwords.

# C.3.1 Adding User Accounts

- Sign in as the administrator.
- Select About > Users in BD Accuri C6 Software.

**NOTE:** The *Users* menu option is only visible to the administrator.

In the Users dialog box, click on Add New User.

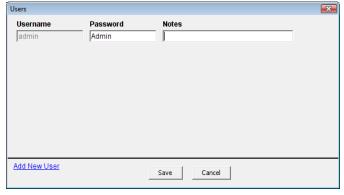


Figure C-4. The Users Information Box

• Type the *Username*, *Password* and *Notes* (optional) in the blank text boxes. Notes are only visible in the Users dialog box.

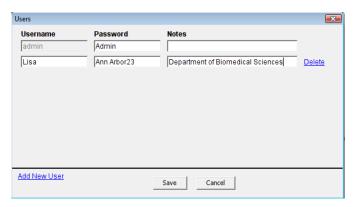


Figure C-5. Adding a New User

Click on the Save button to save changes.

# C.3.2 Deleting User Accounts

- Select About > Users in BD Accuri C6 Software.
- Click on Delete next to the user account to be removed.

**NOTE:** The administrator's account cannot be deleted.



Figure C-6. Deleting a User

· Click on the Save button to save changes.

# C.3.3 Changing a Password

The administrator can change a user's password at any time.

To change a user password:

- Select About > Users in BD Accuri C6 Software.
- Delete the text in the Password field and type a new password.
- Click on the Save button.

# C.4 Signing In and Signing Out

Once tracking is set up, users must sign in for each session.

To sign in to BD Accuri C6 Software:

- Open BD Accuri C6 Software.
- When prompted, type the *Username* in the text box.



Figure C-7. Username and Password Dialog Box

- Type the Password in the text box.
- Click on the OK button
- · Use BD Accuri C6 Software as usual.

BD Accuri C6 Software will continue to log time until the user signs out.

To sign out of the software:

In BD Accuri C6 Software, select File > Quit.

# C.5 Restoring a Forgotten Administrator Password

The password file is encrypted and contains all username and password information created in the User Logging feature. If the Administrator password is forgotten, the password file can be deleted and the Administrator account recreated.

**CAUTION:** This procedure deletes *all* user names and passwords; each account must be manually recreated.

To restore the administrator password:

- If BD Accuri C6 Software is running, shut it down.
- Navigate to the CytometerSupportFiles folder (typically in the root directory of the BD Accuri C6 Software computer).
- · Locate the Password file and delete the file.
- Follow the procedures in sections C.2 and C.3.1.

# **C.6 Monitoring User Activity**

Each time a user signs into, or signs out of, BD Accuri C6 Software, an entry is made in the userUsage Log. The userUsage log is a .csv file.

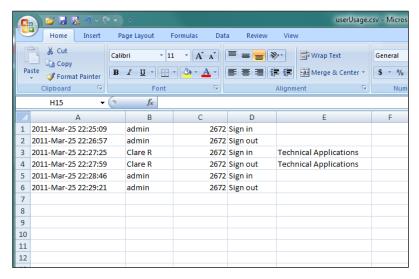


Figure C-8. userUsage Log

The userUsage Log contains the following information:

- Date of a sign in/sign out
- Time of a sign in/sign out
- Username of the operator
- Serial number of the cytometer
- Type of activity (sign in or sign out)

To view the userUsage log:

 Navigate to the CytometerSupportFiles folder (typically in the root directory of the BD Accuri C6 Software computer) and open the folder.

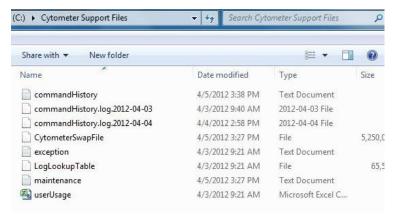


Figure C-9. CytometerSupportFiles Folder

Open the userUsage.log file.

# APPENDIX D BD ACCURI C6 ANALYSIS SOFTWARE

The optional BD Accuri C6 Analysis software allows analysis of BD Accuri C6 Software files on a computer that does not have a connection to the cytometer. BD Accuri C6 Analysis is available in two forms:

- PN 653122—BD Accuri C6 Analysis Software for PC or Mac
- PN 653123—BD CSampler Analysis Software for PC or Mac

BD Accuri C6 Analysis Software has all the functionality of BD Accuri C6 Software, with the exception of instrument control functions. Thresholds and compensation can still be set. BD Accuri C6 Analysis Software is not intended for operating the cytometer.

#### MINIMUM SYSTEM REQUIREMENTS

PC version:

Intel Core Duo processor, 2.8 GHz

2 GB RAM

**CD-ROM Drive** 

Windows® XP (Service Pack 2), Windows Vista, or Windows 7 (recommended)

Language selection: US or UK English

5 GB hard disk space

Display resolution 1280 x 1024 or higher

· Mac version:

Intel or PowerPC processor, 1.66 GHz (Intel processor recommended)

1 GB RAM

**CD-ROM Drive** 

Mac OS X software

Language selection: US or UK English

5 GB hard disk space

Display resolution 1152 x 864 or higher (1280x1024 recommended)

### SOFTWARE OPERATION

See the relevant chapters in this guide for operational information.

# APPENDIX E SELECTABLE LASERS

The detectors and lasers of the standard BD Accuri C6 flow cytometer operate in a predefined configuration: detectors FL1, FL2, and FL3 read blue laser-excited fluorescence emissions and detector FL4 reads red laser-excited emissions. This configuration is referred to as 3 blue 1 red.

The Selectable Lasers Module (PN 653126) allows operation of the cytometer in two alternate configurations which significantly expands the fluorochrome combinations that can be analyzed (see section E.4 for examples):

- 2 blue 2 red:
  - FL1 and FL2 read blue laser-excited emissions FL3 and FL4 read red laser-excited emissions
- 4 blue: All 4 detectors read blue laser-excited emissions

### Components Supplied:

- Selectable Lasers Activation Key (software)
- Three optical filters:
  - 780/60 BP (PN 653187)
  - 610/20 BP (PN 653186)
  - 630/30 BP (only available with Selectable Lasers Module)

**NOTE:** The Selectable Lasers upgrade requires prior installation of BD Accuri C6 Software or BD CSampler Software.

# **E.1** Installing the Selectable Lasers Module

- Verify that BD Accuri C6 Software or BD CSampler Software version 227.4 or above is loaded on the computer. The Selectable Lasers Module can also be installed on any computer where BD Accuri C6 Analysis Software (version 227.4 or higher) is installed.
- Copy the Selectable Lasers Activation Key from the installation CD (or flash drive) to the desktop.
- Double-click on Selectable Lasers Activation Key icon.



Figure E-1. Selectable Lasers Installer

- In the installation wizard, choose the directory in which to install the Activation Key. The location depends on where BD Accuri C6 Software or BD CSampler Software has been installed. In most cases, the location will be in one of the following places:
  - C:\Program Files\BD Accuri\BD Accuri C6 Software\ActivationKeys
  - C:\Program Files\BD Accuri\BD CSampler\ActivationKeys

**NOTE:** If BD Accuri C6 Software or BD CSampler Software are installed on the same computer, install the Selectable Lasers Activation Key twice: once in the BD Accuri C6 Software directory, and again in the BD CSampler Software directory.

Click on the browse button ... to navigate to the correct location.

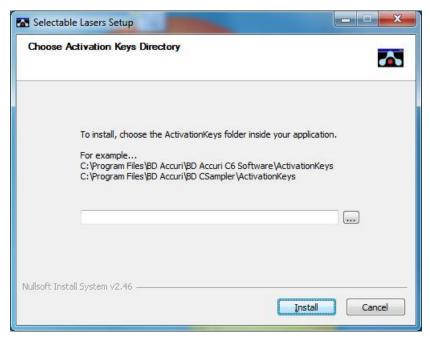


Figure E-2. Install Wizard: Choose Activation Keys Directory

• Click on the Install button.

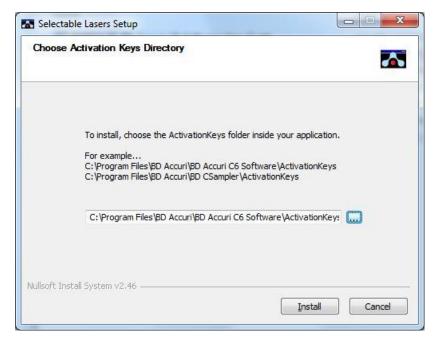


Figure E-3. Install Wizard: Install the Module

• After installation, click on the Close button.

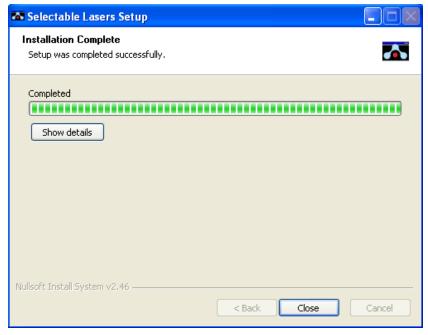


Figure E-4. Install Wizard: Setup Completed Successfully

• Open BD Accuri C6 Software. The Selectable Lasers controls are displayed next to the BD Accuri C6 Software traffic light message.

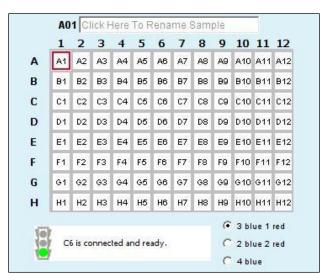


Figure E-5. Selectable Laser Controls in the BD Accuri C6 Software Workspace

 Depending on the laser configuration to be used, the standard optical filters may need to be replaced with one or more of the filters that are supplied with the Selectable Lasers software. See section E.4 for details.

# **E.2** Validating Proper Function After Installation

After installing the Selectable Lasers Module, verify that the module is operating properly. To verify the Selectable Lasers function:

- Leave the standard optical filters in place.
- With BD Accuri C6 Software open and the cytometer powered on, select the 3 blue 1 red laser configuration.



Figure E-6. Selectable Laser Options: 3 Blue 1 Red

- Set the run limit to 30,000 events in all and set the fluidics rate to slow.
- Collect data files for the 8- and 6-Peak Validation Beads (PN 653144 and PN 653145).
- Confirm that the cytometer is operating within specification (see chapter 2).

• Select the 2 blue 2 red option and collect data using the 6-Peak Validation Beads (PN 653145) into a new data well.

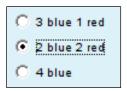


Figure E-7. Selectable Laser Options: 2 Blue 2 Red

 Select the 4 blue option and collect data using the 8-Peak Validation Beads (PN 653144) into a new data well.



Figure E-8. Selectable Laser Options: 4 Blue

Confirm proper operation of the 2 blue 2 red configuration by comparing the 6peak bead distributions for FL3 and FL4 to those obtained with the standard 3
blue 1 red configuration. Results for the 2 blue 2 red selection should show 6
peaks for both FL3 and FL4.

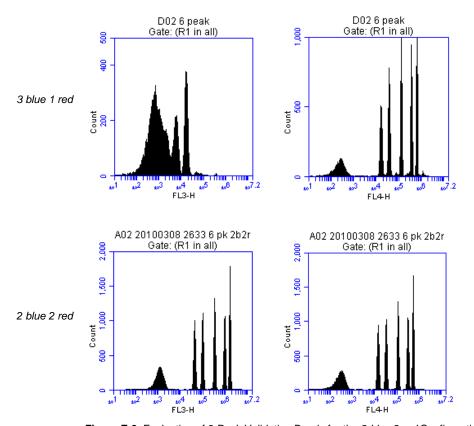


Figure E-9. Evaluation of 6-Peak Validation Beads for the 2 blue 2 red Configuration

 Confirm proper operation of the 4 blue configuration by comparing the 8-peak distributions for FL3 and FL4 to those obtained with the standard 3 blue 1 red configuration. Results for the 4 blue selection should show similar peak profiles for FL3 and FL4.

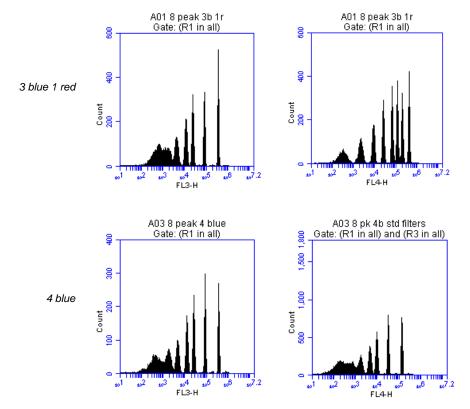


Figure E-10. Evaluation of 8-Peak Validation Beads for the 4 Blue Configuration

# **E.3** Annotating Selected Laser Configuration

The Selectable Lasers Module radio buttons show the *most recently* selected configuration only. BD Accuri C6 Software does not retain this setting for each well. Therefore, it is recommended to annotate the data wells in the naming field of the software to indicate the laser configuration used during data collection of each well, especially when using alternate configurations. For example, name a sample "HPB 4b" to indicate that the *4 blue* option was selected during data collection.

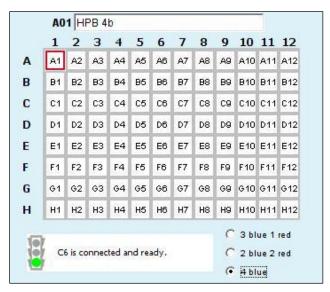


Figure E-11. Renamed Data Well to Indicate 4 Blue Configuration

Laser configurations are *not saved* in the BD Accuri C6 Software data file or with a BD Accuri C6 Software template. Any previously saved BD Accuri C6 Software file will default back to the *3 blue 1 red* option when it is opened.

The last laser configuration used to collect a given well of data is written in the FCS file header using the custom keyword **#LASERCONFIGURATION**. FCS file headers can be viewed by opening an exported FCS file in a text editor (such as Microsoft Notepad).

# **E.4** Optical Filter Placement

Due to the unique optical layout of the cytometer, it is critical that any optional filters used with the Selectable Lasers Module are placed in the proper position for optimal performance.

**WARNING:** The 630/30 bandpass filter provided with the Selectable Lasers Module should only be used when operating in the *4 blue* configuration. Using the 630/30 filter when operating in any other configuration may damage the corresponding detector due to unfiltered red laser signal.

Use the tables below as a guide to optical filter placement for various fluorochrome combinations:

Table E-1. 3 Blue 1 Red: Configuration 1 (Standard Filters)

Detector	Filter	Fluorochrome
FL1	530/30 or 533/30	FITC, GFP, CFSE
FL2	585/40	PE, PI PE-Texas Red
FL3	670 LP	PerCP-Cy5.5 PE-Cy5, PE-Cy7

FL4	675/25	APC, Alexa-647 PE-Cy5
-----	--------	--------------------------

Table E-2. 3 Blue 1 Red: Configuration 2

Detector	Filter	Fluorochrome
FL1	530/30 or 533/30	FITC, GFP, CFSE
FL2	585/40	PE
FL3	610/20	PI PE-Texas Red
FL4	675/25	APC, Alexa-647 PE-Cy5

**NOTE:** When operating in the *3 blue 1 red* configuration, place either the 610/20 bandpass or the 670 LP filters in position FL3. For best results when analyzing PE and PE-Texas Red (PE-TR) simultaneously, select the filters with the following signal-intensity considerations in mind:

- PE-bright, PE-TR-moderate to bright: FL3 = 670 LP
- PE-dim to moderate, PE-TR any level: FL3 = 610/20
- PE-bright, PE-TR dim: may be difficult to separate; consider using the *4 blue* configuration with a 630/30 in FL4 to detect PE-TR.

Table E-3. 2 Blue 2 Red Configuration

Detector	Filter	Fluorochrome
FL1	530/30 or 533/30	FITC, GFP, CFSE
FL2	585/40	PE, PI PE-Texas Red
FL3	780/60	APC-Cy™7 (and equivalents)
FL4	675/25	APC (and equivalents)

Table E-4. 4 Blue: Configuration 1

Detector	Filter	Fluorochrome
FL1	530/30 or 533/30	FITC, GFP, CFSE
FL2	585/40	PE, PI PE-Texas Red
FL3	780/60	PE-Cy7 (and equivalents)
FL4	675/25	PerCP-Cy5.5 PE-Cy5

Table E-5. 4 Blue: Configuration 2

Detector	Filter	Fluorochrome
FL1	530/30 or 533/30	FITC, GFP, CFSE
FL2	585/40	PE
FL3	780/60	PE-Cy7 (and equivalents)
FL4	610/20	PI
	630/30	PE-Texas Red

Table E-6. 4 Blue: Configuration 3

Detector	Filter	Fluorochrome
FL1	530/30 or 533/30	FITC, GFP, CFSE
FL2	585/40	PE
FL3	675/25	PerCP-Cy5.5 PE-Cy5
FL4	610/20	PI
	630/30	PE-Texas Red

# **E.5 Selectable Laser Application Examples**

# E.5.1 2 Blue 2 Red Configuration Examples

Table E-7. Detector and Filter Configuration for 2 Blue 2 Red, Examples 1 and 2

Detector	Filter	Fluorochrome
FL1	530/30 or 533/30	FITC
FL2	585/40	PE
FL3	780/60	APC-Cy7
FL4	675/25	APC

## Example 1: 2 Blue 2 Red for Simultaneous FITC, PE, APC and APC-Cy7

The following images illustrate a 4-color analysis of CD3<sup>+</sup>CD4<sup>+</sup> cells. Human peripheral blood (HPB) was stained with CD3-APC-Cy7, CD4-APC, CD45RA-FITC, and CD45RO-PE.

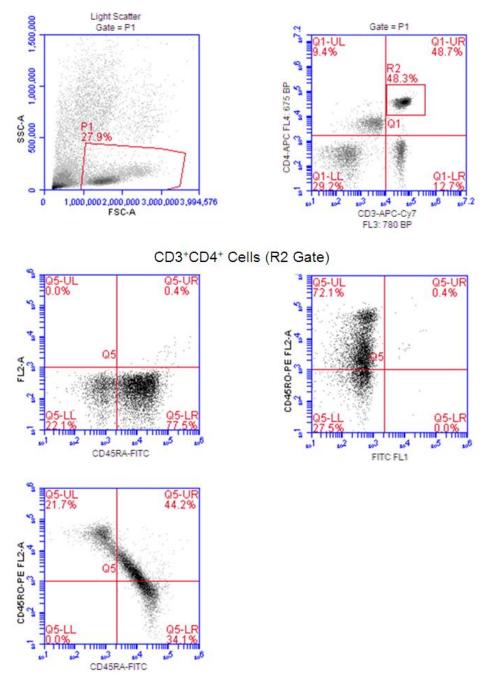


Figure E-12. Gating Example Using the 2 Blue 2 Red Configuration

# Example 2: 2 Blue 2 Red with BD Cytometric Cytokine Bead Array (CBA) 30 Plex Bead Mixture (BD 51-9004679)

The following images show the results for bead separation using the entire 30 plex bead mixture

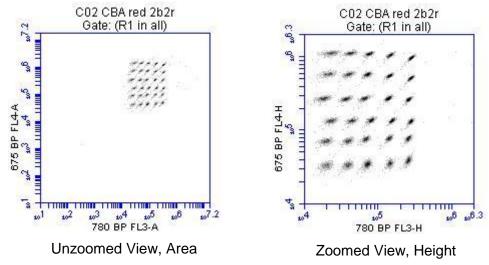


Figure E-13. BD Accuri C6 Software (PN 653126).

# E.5.2 4 Blue Configuration Examples

## Example 1: 4 Blue FITC, PE, PE-Cy5, PE-Cy7: Human Peripheral Blood Subsets

Table E-8. Detector and Filter Configuration for Selectable Lasers 4 Blue Example 1

Detector	Filter	Fluorochrome
FL1	530/30 or 533/30	FITC
FL2	585/40	PE
FL3	780/60	PE-Cy7
FL4	675/25	PE-Cy5

The following images show CD45-FITC vs. SSC gating of lymphocytes (zoomed in, left; zoomed out, right). HPB was stained with CD45-FITC, CD8-PE, CD3-PE-Cy5, and CD4-PE-Cy7).

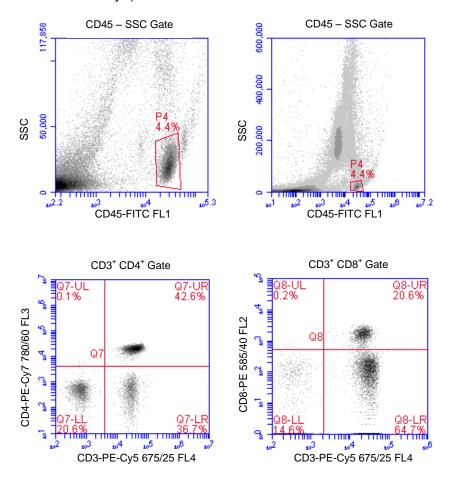


Figure E-14. 4 Blue Configuration and the Optional 780/60 Optical Filter at Detector FL3

# Example 2: 4 Blue FITC, PE, PE-Cy5, PE-Texas Red: Human Peripheral Blood Subsets

Table E-9. Detector and Filter Configuration for Selectable Lasers 4 Blue, Example 2

Detector	Filter	Fluorochrome
FL1	530/30 or 533/30	FITC
FL2	585/40	PE
FL3	675/25	PE-Cy5
FL4	630/30	PE-Texas Red

The following images show the results for example 2. The color compensation value to correct the spillover of PE (FL2) into the PE-Texas Red (FL4) detector may be in the range of 70% to 90%.

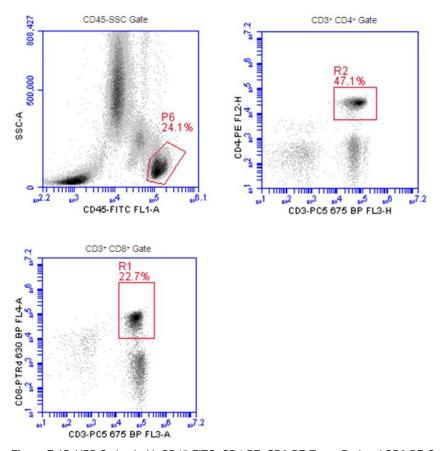


Figure E-15. HPB Stained with CD45-FITC, CD4-PE, CD8-PE-Texas Red and CD3-PE-Cy5

# APPENDIX F ENHANCED ANALYSIS FEATURES

The Enhanced Analysis module includes several advanced functions in BD Accuri C6 Software:

- · Live (or "dump") gating
- Plot and region renaming
- Event coloring
- Vector scalable graphics
- Batch analysis

**NOTE:** Enhanced Analysis is compatible with BD Accuri C6 Software version 264 and higher.

# F.1 Creating a Live Gate

Live gating, also called dump gating, is used to exclude events outside a designated region from being displayed and stored as part of the c6 or fcs file. Any region may be used as a live gate. This selection can be remembered as part of the template if desired.

**CAUTION:** Once a live gate is executed, there is no way of recapturing any excluded data.

To create a live gate:

- Create the region to use in the live gate in the Collect Tab.
- Click on the *Do not collect events outside* check box and select the region from the drop-down list located under *Run Settings* in the *Collect Tab*.

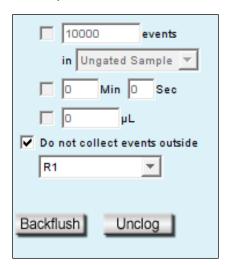


Figure F-1. Creating a Live Gate

# F.2 Renaming Plots and Regions

Plots or regions in a plot can be renamed. In each case the method is the same.

To rename a plot or region:

· Double-click on the plot or region name.

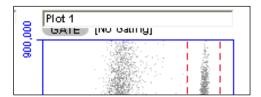


Figure F-2. Enable Renaming of the Plot

In the text field, type in a new name.

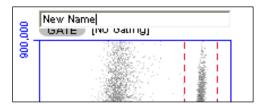


Figure F-3. Type a New Plot Name

· Press Enter.

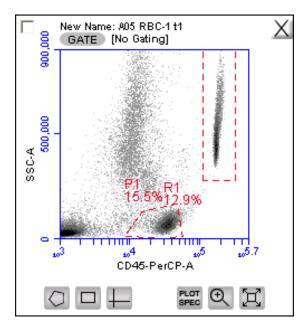


Figure F-4. New Plot Name

# F.3 Coloring Events in a Region

All events within one or more regions can be designated to appear as a specified color within other histogram or dot plots. The event coloring will remain and automatically update during data acquisition.

NOTE: Color will not appear in a density plot.

To color events:

- Create a region in a histogram, dot or density plot.
- Double-click on the region name and click on the small white square to display a color palette. The most commonly used colors appear on the top row.

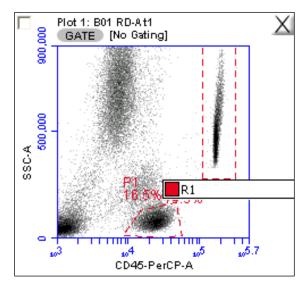


Figure F-5. Select a Region to Color

Select a color.

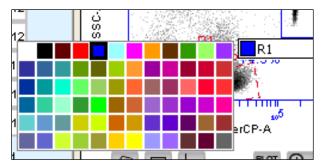


Figure F-6. Select a Color for a Region

 Click outside the region to make the change take effect. Events that fall within the region in other dot and histogram plots are now colored and displayed in the same color.

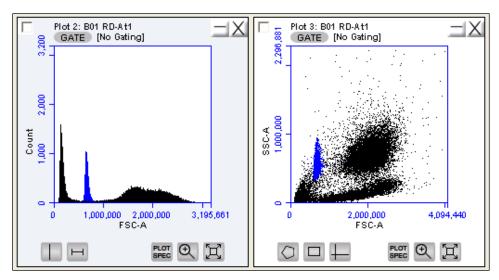


Figure F-7. Colored Events with a Region

# F.4 Creating Publication-Quality Images of Plots

High resolution (publication quality) images are stored in eps format, which is vector scalable.

To create a high resolution image of a plot:

- Select File > Set Plot Drag and Drop Format.
- Select the .eps option in the dialog box.
- Click on the OK button.
- In the BD Accuri C6 Software Workspace, click on a plot and drag it onto the desktop.
- Open an image editing application (e.g., Photoshop) and ensure the image resolution is set to 300 dpi, or higher if required, and import the eps image.
- Save the image.

# F.5 Analyzing Batches of Samples

The *Batch Analysis* tab allows an automated analysis to be performed on multiple samples at the same time. The plot types created during acquisition will appear at the top of the screen. Statistics can be displayed for individual files or all files selected. After analysis, regions can be moved or resized on plots for individual samples without affecting the plots for other samples.

# F.5.1 Viewing the Batch Analysis Tab

The Batch Analysis tab is organized into two major sections:

 Setup panel—Panel on the top half of the window that contains controls for selecting samples and plots for analysis.

 Data display—Large area on the bottom half of the window that shows sample data in plots and in a Statistics Table.

When the *Batch Analysis* tab is opened for the first time, the workspace is empty. To set up the *Batch Analysis* tab, plots can be created or copied from the *Collect or Analyze* tabs.

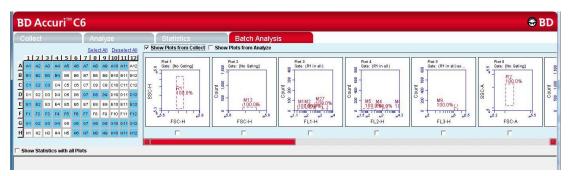


Figure F-8. Batch Analysis Tab Workspace

The following table describes each of the controls and indicators in the *Batch Analysis* tab:

Table F-9. Batch Analysis Tab Controls

Control	Description						
Sample Grid	Matrix laid out in the configuration of a 96-well plate for selecting the samples to analyze.						
	The wells are color-coded:						
	White—Does not contain data.						
	Blue—Contains data.						
	Black check mark—Currently selected for batch analysis.						
Available Plots	Displays plots (without data) that can be included in the batch analysis. Available plots include plots copied from the <i>Collect</i> tab or created in the <i>Analyze</i> tab. See section F.5.2 for details.						
Analysis Pane	Rows of analysis data in table and plot format, based on the plots that were selected for use in analysis.						
	Statistics Tables can be copied into most Microsoft Office compatible programs.						

# F.5.2 Running a Batch Analysis

- · Run a number of samples.
- After running the samples, click on the Batch Analysis tab.
- Select the samples to analyze in the Sample Grid by clicking on the wells.
- Do one or both of the following:

- Enable the Show Plots from Collect check box to make all of the plots from the Collect tab available for batch analysis.
- Enable the Show Plots from Analyze check box to make all of the plots from the Analyze tab available for batch analysis.
- Select the specific plots to include in the batch analysis by enabling the check boxes under those plots.
- To show the statistics associated with samples, do one of the following:
  - Enable the Show Statistics with All Plots check box to show statistics in every row.
  - Enable the *Show Statistics* check box in individual rows to view statistics for specific samples. Regions can be moved and resized (see section 3.7).

## F.5.3 Exporting Data

Plots and statistics can be exported in two ways, either as a PowerPoint file or as an Excel spreadsheet. To export data:

- Do one of the following in the lower right-hand corner of the *Batch Analysis* tab:
  - Click on the Export to PowerPoint button.
  - Click on the Export Stats to Excel button.



Figure F-10. Export Buttons in the Batch Analysis Tab

In the Save dialog box, navigate to the destination folder and type a file name. BD Accuri C6 Software exports the file to the specified location.

# APPENDIX G FCS KEYWORDS

The following tables list the keywords used in BD Accuri C6 Software FCS files.

## STANDARD FCS TAGS IN FCS FILES EXPORTED BY BD ACCURI C6 SOFTWARE

Table G-1. Standard FCS Tags

Table G-1. Standard FCS Tage	S	
FCS Tag	Description	BD Accuri C6 Software
\$FIL	Filename including .fcs extension.	Save file dialog.
\$SMNO	Sample name.	Sample name field above the Well Grid.
\$DATATYPE	The data type of the actual values for each event. It is always "I" for unsigned binary integers.	Unchangeable.
\$MODE	The mode of the data. It is always "L" for list mode where the data is in the order described by the \$Pn keywords.	Unchangeable.
\$BYTEORD	Order in which data bytes are written, least to most significant. It is always 4,3,2,1.	Unchangeable.
\$NEXTDATA	The byte offset for an additional dataset in the file. BD Accuri files always specify 0 since the files only contain 1 dataset.	Unchangeable.
\$PAR	Total number of parameters stored in the dataset. All datasets have 14 parameters.	Unchangeable.
\$PnB	For parameter N, the number of bits for each binary value. The number is always 32 since integers are stored in 32 bits in Java.	Unchangeable.
\$PnR	The range of parameter N. The range for all parameters is always 16777216.	Unchangeable.
\$PnN	The name of parameter N. Parameters are the default values from BD Accuri C6 Software.	Unchangeable.
\$PnE	For parameter N, this denotes if linear or logarithmic amplifiers are used. It is always 0,0 because linear values for data are always saved. This is an optional tag.	Unchangeable.
\$PnS	The name of the fluorescence stain or probe used for parameter N. This tag is used for the custom parameter name.	Rename parameter dialog.
\$ТОТ	Number of objects stored in the data list. The cumulative event total for the sample.	Determined by the amount of events acquired by user.
\$DATE	The date the represented sample was last acquired into DD-MMM-YYYY.	Defined by the beginning time of the most recent acquisition. Set by the computer's clock.

FCS Tag	Description	BD Accuri C6 Software
\$CYT	The name of the cytometer used for the measurement. It is always BD Accuri C6.	Unchangeable.
\$SPILLOVER	The standard tag for color compensation.	Determined by values entered into Color Compensation dialog.
\$TIMESTEP	Hard-coded value of 0.1, which is in seconds.	Unchangeable.
\$PROJ	Represents the name of the workspace, which is also the name of the .c6 file (before the .c6 extension). If exporting an FCS file, this value is the name of the file before the .fcs extension.	Saving as a different file, importing an FCS file, or exporting an FCS file would change this value.
\$BTIM	The beginning time of acquisition of the first event.	Unchangeable.
\$ETIM	The end time of acquisition of the last event.	Unchangeable.
\$VOL	Total volume in nanoliters.	Collecting more events would change the total volume.
\$BEGINSTEXT	Default FCS 3.0 tag to mark the beginning of the text section.	Unchangeable.
\$ENDSTEXT	Default FCS 3.0 tag to mark the end of the text section.	Unchangeable.
\$BEGINANALYSIS	Default FCS 3.0 tag to mark the beginning of the analysis section.	Unchangeable.
\$ENDANALYSIS	Default FCS 3.0 tag to mark the end of the analysis section.	Unchangeable.
\$BEGINDATA	Default FCS 3.0 tag to mark the beginning of the data section.	Unchangeable.
\$ENDDATA	Default FCS 3.0 tag to mark the end of the data section.	Unchangeable.

## **CUSTOM TAGS IN FCS FILES EXPORTED BY BD ACCURI C6 SOFTWARE**

Table G-2. Custom FCS Tags

Table G-2. Custom FCS Tags		
Tag	Description	BD Accuri C6 Software
#BDACCURIDECADESN	The number of decades for parameter N. It is always 7.224719870049579.	Unchangeable.
#BDACCURI4COLORCOMP	A list of the percent values as entered into the color compensation window.	Determined by values entered into Color Compensation dialog.
#BDACCURICAPTUREDDATE	The date of the last time the represented sample was acquired into expressed in milliseconds since Jan 1, 1970.	Defined by the beginning time of the most recent acquisition. Set by the computer's clock.
#PNVIRTUALGAIN	The VirtualGain set for parameter N, where 1.0 means no VirtualGain.	Set by user in the VirtualGain window.
#SAMPLE	Value is either the well code or the sample rename (if one exists).	Can be changed by renaming the sample in the application. If no rename exists, importing an FCS file into a different well than the one collected in will change this value.
#ATIM	Cumulative acquisition time.	Can be changed by acquiring more data.
#SPACERS	Used to pad the text, analysis, and data sections of the FCS file.	Unchangeable.
#LASERCONFIGURATION	Designates selected mode of operation when FCS file was collected: 3 blue 1 red, 2 blue 2 red, or 4 blue.	Changes if a different mode is selected and then data is added to a well.
#PnMaxUsefulChannelDNA	The data value for parameter N that is the 95 <sup>th</sup> percentile.	Defined by data values.

## FCS 3.0 DEFINED TAGS NOT IN BD ACCURI C6 SOFTWARE EXPORTED FCS FILES

Table G-3. FCS Tags Not in BD Accuri C6 Software Exported Files

Tag	Description
\$COMP	Amount of fluorescence compensation employed during collection. This replaces the \$DFCiTOj tag from FCS 2.0.
\$TIMESTEP	Absolute measure of time used in kinetic analysis.
\$UNICODE	Enables usage of certain keywords in non-English languages. This is optional.
\$PnE	For parameter N, this denotes if linear or logarithmic amplifiers are used. BD Accuri C6 Software always uses 0,0 because BD Accuri C6 Software always saves linear values for the data. Mandatory in 3.0.
\$PnN	When time is collected, the keyword value of the time- parameter name must now be the string "TIME".
CRC	An optional 16-bit Cyclic Redundancy Check has been added to the end of each dataset.

# APPENDIX H EXAMPLE .CSV FILE

The following table is an exported bead log file in .csv format.

Each row represents the parameter data associated with a single event. The number of rows will equal the total number of events in the data file. A data well that contains 125,000 events also contains 125,000 rows of data in the spreadsheet display of the .csv data. The Time value is rounded to the nearest second during data export. The Width parameter is that for the primary threshold parameter.

Table H-1. Example .csv File

	Α	В	С	D	E	F	G	Н	T.	J	K	L	M	N
1	FSC-A	SSC-A	FL1-A	FL2-A	FL3-A	FL4-A	FSC-H	SSC-H	FL1-H	FL2-H	FL3-H	FL4-H	Width	Time
2	747359	206049	785	147	0	423	1883688	704036	486	60	161	382	82	137
3	308831	36583	390	221	456	0	787489	127133	274	114	438	383	75	137
4	656756	159245	685	179	870	286	1935215	555473	443	114	780	652	78	137
5	519654	138403	746	329	225	763	1175444	463213	501	152	118	506	86	137
6	818135	257287	767	72	1068	691	2171916	876805	591	0	586	629	80	137
7	257396	19821	203	254	75	76	688493	71597	110	96	182	298	73	137
8	268765	16276	168	5	0	26	742169	58897	46	55	275	362	71	137
9	741749	183316	329	145	517	526	2209348	664629	200	48	431	458	77	137
10	74788	19503	295	204	70	194	279103	76864	154	70	0	278	44	137
11	310347	28687	405	213	513	206	811460	101016	259	79	345	631	75	137
12	676065	201903	585	500	1374	775	1755862	709407	432	208	702	594	81	137
13	611517	252169	788	242	400	0	1461590	889724	427	177	274	229	82	137
14	321267	42064	149	194	244	660	823497	146225	319	46	314	732	76	137
15	301525	56315	314	133	442	118	734258	189713	150	56	414	442	78	137
16	787700	328238	926	92	1402	15	2011475	1183340	798	66	823	454	82	137
17	712542	188615	567	84	376	0	1939139	664650	415	48	343	234	79	137
18	940993	240525	871	224	122	0	2740219	850548	669	85	121	252	79	137
19	93345	18538	65	110	786	378	340737	73594	34	14	668	293	49	137
20	1120179	898131	6328	1672	1878	3020	3199136	3339490	5642	1524	1869	2551	80	137
21	735806	205623	342	158	776	565	2127351	755878	229	131	364	480	76	137
22	93556	28907	425	178	1106	217	312696	108595	204	66	464	318	51	138
23	539263	150426	657	220	626	435	1257974	519183	631	124	330	390	82	138
24	390039	401617	720	240	1270	542	1519975	1555340	604	108	747	613	57	138
25	656006	159032	570	198	433	0	1809412	567177	340	205	243	380	79	138
26	749099	186852	526	177	437	374	2095595	670857	287	81	187	407	79	138
27	249477	24622	598	267	757	0	640125	90865	424	97	486	378	74	138

# APPENDIX I TECHNICAL SUPPORT

# For Technical Support, contact

#### USA

Accuri Cytometers, Inc. 173 Parkland Plaza Ann Arbor, MI 48103 USA Phone: +1.734.994.8000 Fax: +1.734.994.8002

Email: <u>CustomerService@AccuriCytometers.com</u> Email: <u>TechSupport@AccuriCytometers.com</u>

## • Europe, Middle East and Africa

BD Biosciences Erembodegem-Dorp 86 9320 Erembodegem Belgium

Phone: +32 53 720 882

Customer Support: <a href="mailto:bd\_accuri@europe.bd.com">bd\_accuri@europe.bd.com</a>
Technical Support: <a href="mailto:techsupport@europe.bd.com">techsupport@europe.bd.com</a>

## Canada

BD Biosciences 2100 Derry Rd West, Suite 100 Mississauga, Ontario Canada L5N 0B3

Email: canada@bd.com

## **Direct Phone Numbers**

Austria: 0810 101 807

Belgium: 078 166 050

Canada: 1.800.268.5430

Denmark: 8025 0622

Finland: 0800 915582

France: 0811 290 069

Germany: 0180 100 1732

**Ireland: 1850 930 396** Italy: 840 999 926

Netherlands: 0900 0400 142

Norway: 800 17 382 **Poland: 00800 121 4744** 

#### Regional Offices bdbiosciences.com/offices

Asia Pacific Singapore Tel 65.6861.0633

Fax 65.6860.1590 **China** 

Tel 86.21.3210.4610 Fax 86.21.5292.5191 India

Tel 91.124.2383566 Fax 91.124.2383224/25/26 Australia/New Zealand Australia

Toll Free: 1800 656 100
Tel 61.2.8875.7000
Fax 61.2.8875.7200
bd\_anz@bd.com
New Zealand

Toll Free: 088 572.468 Tel 64.9.574.2468 Fax 64.9.571.2469 bd\_anz@bd.com Europe BD Biosciences

Tel 32.2.400.98.95 Fax 32.2.401.70.94 bdbiosciences.com/eu help.biosciences@europe.bd.com

Canada

BD Biosciences
Toll Free 888.259.0187
Tel 905.542.8028
Fax 888.229.9918

bdbiosciences.com/ca canada@bd.com Japan

Nippon Becton Dickinson Toll Free 0120.8555.90

Tel 81.24.593.5405 Fax 81.24.593.5761 United States BD Biosciences

Toll Free 877.232.8995 Fax 800.325.9637 bdbiosciences.com answers@bd.com

For Research Use Only. Not for use in diagnostic or therapeutic procedures. Class 1 Laser Product.

BD Accuri, BD Accuri C6 Software, BD Accuri CSampler Software, BD CSampler, VirtualGain, BD Accuri C6 Flow Cytometer and "Science is hard. Flow cytometry should be easy" are trademarks of Becton Dickinson and Company. ©2012