
Operating Instructions
Bedienungsanleitung
Mode d'emploi
Manual de Instrucciones
Istruzioni d'uso

METTLER TOLEDO

METTLER TOLEDO
RS-P42 Printer
Drucker RS-P42
Imprimante RS-P42
Impresora RS-P42
Stampante RS-P42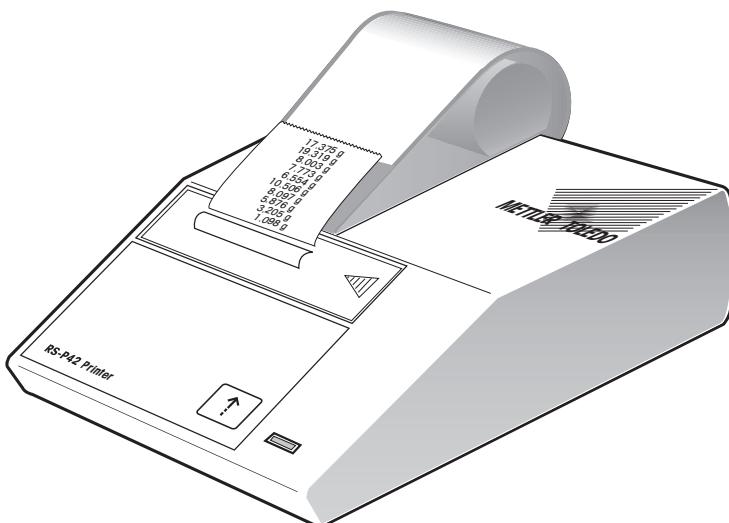

Page 3

English

Seite 10

Deutsch

Page 17

Français

Página 24

Español

Página 31

Italiano

Content

1	Introduction	4
1.1	What are the capabilities of the RS-P42 Printer	4
2	Startup procedure	4
2.1	Safety measures	4
2.2	Adjust Baudrate	5
2.3	Inserting paper	5
2.4	Attaching the printer	5
3	Operation	6
4	Printer test	6
5	Maintenance	6
5.1	Changing paper	6
5.2	Replacing the ribbon	7
5.3	Cleaning	7
6	Malfunctions	7
7	Consumables	8
8	Technical data	8

1 Introduction

1.1 What are the capabilities of the RS-P42 Printer

The RS-P42 is a rugged and simple to operate dot matrix printer for METTLER TOLEDO balances with an RS232C interface. The printout on normal paper fulfills the requirements of modern quality assurance systems (GLP, GMP, ISO 9001, etc.). These define, among other things, the principles and duties for the user of measuring instruments or inspection, measuring and test equipment. The RS-P42 complies with the principle of traceability of measurements by precisely documenting the origin of data, as well as calibration procedures. It fulfills the obligation to archive measurement results by its thermally stable and lightfast print. On connection to METTLER TOLEDO balances, the RS-P42 prints all data sent by the balance without restriction, that is not only weight values and calibration procedures, but also the results of all applications, built into the balance.

2 Startup procedure


2.1 Safety measures

The RS-P42 has been tested for the connections and intended purposes documented in these Operating Instructions. However, this does not absolve you from the responsibility of performing your own tests of the product supplied by us regarding its suitability for the methods and purposes you intend to use it for. You should therefore observe the following safety measures.


- The RS-P42 Printer may be used only indoors and on no account in hazardous areas.
- The electronics of the RS-P42 Printer are protected against the ingress of splashing water and dust. However, as the printer housing is not tight the printer should not be used in the presence of liquids.
- Comply with cleaning instructions (see item 5.3), protect power plug against wet conditions.
- Never open the printer housing. Should you ever have problems with your printer, please contact your responsible METTLER TOLEDO dealer.

2.2 Adjust Baudrate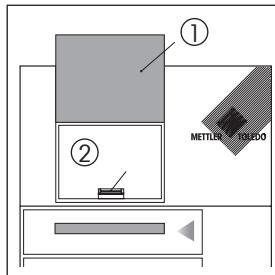

- Open cover of paper compartment (1) and remove paper roll.
- Press plastic tab (2) backward and open the compartment cover.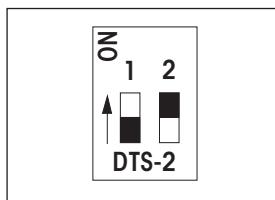

Baudrate


1200/8/N/None
2400/7/E/None

Dip-Switch

1	2
on	off
off	off (Factory setting)

Important: To ensure that special characters (e.g. "°C") are printed out correctly on METTLER TOLEDO printers, the **balance and printer** must be set to **1200/8/N/None**.


2.3 Inserting paper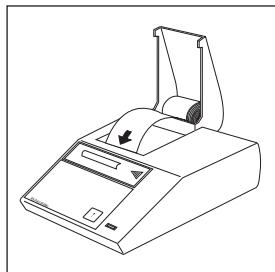

- Feed paper through the slot in the printing unit in the direction of the arrow.
- Switch on printer with toggle switch at the rear.
- Press and hold until sufficient paper has been fed in.

English

2.4 Attaching the printer

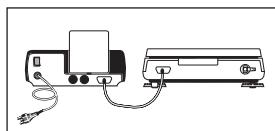- Attach connection cable of the RS-P42 to the printer interface of the instrument and to the printer.
After configuring the attached instrument for the RS-P42, the printer is ready for operation.

3 Operation

- Pressing the printer key advances the paper 1 line (line feed).
- Pressing the balance key with the symbol prints out the stable weight value.
As long as the balance does not show a stable weight value, nothing will be printed out.

4 Printer test

The RS-P42 is equipped with a self-test which automatically prints out the character set of the printer.

Starting the test

- Switch on RS-P42 and at the same time press and hold for a few seconds.

Terminating the test

- Switch off the RS-P42.


5 Maintenance

Under normal ambient conditions, the RS-P42 Printer requires practically no maintenance thanks to its rugged construction. The METTLER TOLEDO service department is available should servicing be required. Please ask your responsible METTLER TOLEDO dealer for details.


5.1 Changing paper

Insertion of a new roll of paper is described at item 2.3. Order number for paper, see item 7.

5.2 Replacing the ribbon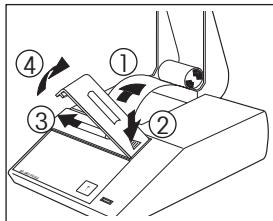

- Pull paper out of printing unit (1).
- Remove ribbon cover (2, 3, 4).
- To remove ribbon cartridge, press in direction of arrow (5).
- Insert new ribbon and if need be, take up slack with tensioning wheel (6).
- Refit ribbon cover.
- Insert paper and feed through slit in the printing unit as far as it will go.
- Press and hold until sufficient paper has been fed in.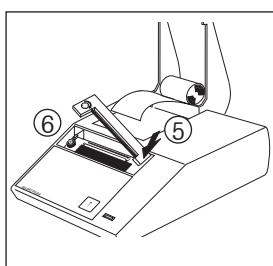

5.3 Cleaning

As the printer housing is made of top grade, resistant materials, all commercially available cleaning agents may be used.

English

6 Malfunctions

Malfunction / Message	Possible cause	Rectification
Printout unreadable	<ul style="list-style-type: none"> • Ribbon worn / entangled • Service life of printing unit at an end 	<ul style="list-style-type: none"> • Change / tension ribbon • Please contact METTLER TOLEDO service
Green pilot lamp does not light up	<ul style="list-style-type: none"> • Printer on • No power, fuse defective 	<ul style="list-style-type: none"> • Please contact METTLER TOLEDO service

7 Consumables

	Type / Order number	Standard
Paper roll, set of 5	72456	2 piece
Paper roll, adhesive, set of 3	11600388	1 piece
Ribbon cartridge, black, minimum order quantity: 2	65975	1 piece
Connection cable for: METTLER TOLEDO balances with an RS232C interface	11101051	1 piece
Spare printing unit (may be changed only by METTLER TOLEDO service!)	600133	–
Power line fuse 230 V (T100L250V) or 115 V (T200L250V) (may be changed only by METTLER TOLEDO service!)		commercially available

8 Technical data

Printer


Printing unit	Dot matrix printer 5x9 dots, 24-character line length,
Epson	IBM character table No. 4
Printing speed	1 line per second
Ribbon cartridge	Exchangeable, black
Paper roll	Normal paper, 58 x dia. 51 mm, integrated in housing, commercial size

Printer functions

Printing, lightfast and thermally stable (GLP, GMP,
ISO 9001), line feed, self-test

RS232C-Interface

The RS-P42 has an RS232C interface for the attachment of METTLER TOLEDO instruments and peripherals.

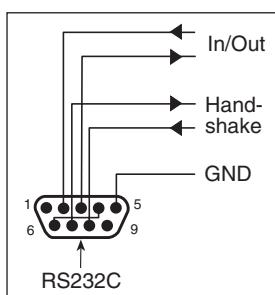- 9-pin male connector (pin assignment as IBM laptop)
- Matching to other device (transmission parameters) see item 2.2
- The hardware handshake lines are active when "None" has been selected in switch position HANDSHAKE

Designed test conditions

Power line voltage, frequency	115 V or 230 V, admissible fluctuations –20% +15%, 50/60 Hz
Power consumption	140 mA or 70 mA
Altitude	4000 m
Temperature range	0...40 °C
Humidity	Max. relative humidity 80% for temperatures to max. 31 °C, linear decreasing to 50% at 40 °C
Installation category	II
Pollution degree	2
Dimensions / Weight	W x D x H = 157 x 210 x 85 mm / net 1.5 kg (incl. paper roll)

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To protect your METTLER TOLEDO product's future:
METTLER TOLEDO Service assures the quality, measuring accuracy and preservation of value
of all METTLER TOLEDO products for years to come.
Please send for full details about our attractive terms of service.
Thank you.

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P11780561

Subject to technical changes and to the availability
of the accessories supplied with the instruments.