

Operating Manual

Translation of the original operating manual

KB / KB-UL (E4), KB / KB-UL (E6) Cooling Incubators

with compressor technology and program control

| Model | Model version | Art. No. |
|----------------|---------------|----------------------|
| KB 53 (E4) | KB053-230V | 9020-0199, 9120-0199 |
| KB 53-UL (E4) | KB053UL-120V | 9020-0302, 9120-0302 |
| KB 115 (E4) | KB115-230V | 9020-0397, 9120-0397 |
| KB 115-UL (E4) | KB115UL-120V | 9020-0398, 9120-0398 |
| KB 240 (E6) | KB240-230V | 9020-0202, 9120-0202 |
| KB 240-UL (E6) | KB240UL-120V | 9020-0304, 9120-0304 |
| KB 400 (E6) | KB400-230V | 9020-0203, 9120-0203 |
| KB 400-UL (E6) | KB400UL-120V | 9020-0305, 9120-0305 |
| KB 720 (E6) | KB720-230V | 9020-0204, 9120-0204 |
| KB 720-UL (E6) | KB720UL-240V | 9020-0306, 9120-0306 |

BINDER GmbH

- ► Address: Post office box 102, 78502 Tuttlingen, Germany ► Phone: +49 7462 2005 0
- ► Fax: +49 7462 2005 100 ► Internet: http://www.binder-world.com
- ► E-mail: info@binder-world.com ► Service Hotline: +49 7462 2005 555
- ► Service Fax: +49 7462 2005 93 555 ► Service E-Mail: customerservice@binder-world.com
- ► Service Hotline USA: +1 866 885 9794 or +1 631 224 4340 x3
- ► Service Hotline Asia Pacific: +852 390 705 04 or +852 390 705 03
- ► Service Hotline Russia and CIS: +7 495 988 15 16

Contents

| 1. | SAFETY | 6 |
|--------------|--|-----|
| 1.1 | Personnel Qualification | |
| 1.2 1.3 | Operating manual Legal considerations | |
| 1.4 | Structure of the safety instructions | |
| 1.4 | I.1 Signal word panel | 7 |
| 1.4 1.4 | | |
| 1.4 | | |
| 1.5 | Localization / position of safety labels on the chamber | 9 |
| 1.6 1.7 | Type plate General safety instructions on installing and operating the chambers | |
| 1.8 | Intended use | |
| 1.9 | Foreseeable Misuse | .14 |
| 1.10 1.11 | Residual Risks Operating instructions | |
| 1.12 | Measures to prevent accidents | |
| 2. | CHAMBER DESCRIPTION | 18 |
| | | |
| 2.1 2.2 | Chamber overview Instrument panel | |
| 2.3 | Lateral control panel with main power switch and optional equipment – KB / KB-UL 240 / 400 / 7 | |
| 0.4 | (E6) | |
| 2.4 | Chamber rear – KB / KB-UL 53 / 115 (E4) | .22 |
| 3. | COMPLETENESS OF DELIVERY, TRANSPORTATION, STORAGE, AND INSTALLATION | 23 |
| 3.1 | Unpacking, and checking equipment and completeness of delivery | |
| 3.2 | Guidelines for safe lifting and transportation | |
| 3.3 3.4 | Storage Location of installation and ambient conditions | |
| 4. | INSTALLATION OF THE CHAMBER | |
| 4.1 | Spacers for wall distance – KB / KB-UL 240 / 400 / 720 (E6) | 27 |
| 4.2 | Mounting the flexible tilt protection kit | |
| 4.3 | Electrical connection | .29 |
| 5. | START UP | 30 |
| 6. | FUNCTIONAL OVERVIEW OF THE T4.12 CHAMBER CONTROLLER | 30 |
| 6.1 | Menu structure | .31 |
| 6.1 | | |
| 6.1 6.1 | | |
| 6.2 | Operating modes | |
| 6.2 | Activating the "control off" mode or change to "fixed value" operating mode | .34 |
| 6.3 6.4 | Performance during and after power failure Information | |
| 7. | CONFIGURATION OF OPTIONAL EQUIPMENT | |
| 7.1 | Turning on / off the optional interior socket | |
| 7.2 | Switching on or off the optional zero-voltage relay control outputs | |
| 7.3 | Functional test of the optional alarm output | .39 |
| 7.4 | Switching on or off the optional object temperature display | .39 |

| 8. | SET-POINT ENTRY IN "FIXED VALUE" OPERATING MODE | 40 |
|--------------|--|----|
| 8.1 | Set-point ranges | 40 |
| 8.2 | Entering the set-points via "quick menu" | |
| 8.3 | Entering the set-points via general menu | |
| 9. | TIME PROGRAMS | 43 |
| 9.1 | Starting and running an existing time program | 45 |
| 9.2 | Cancelling a running time program | |
| 9.3 | Creating a new time program | |
| | 3.1 Section handling 3.2 Temperature setpoint | |
| | 3.3 Section duration | |
| | 3.4 Repeating one or several sections within a time program | |
| 9.3 | 3.5 Tolerance range | 54 |
| | 3.6 Set-point ramp and set-point step modes | |
| | 3.7 Switching on or off the optional zero-voltage relay outputs | |
| | Calling up the next parameter Saving the time program and leaving the program editor | |
| 9.4 | Program interruption | |
| 9.5 | Deleting a time program | |
| 10. | WEEK PROGRAMS | 64 |
| - | | - |
| 10.1 | Starting and running an existing week program | |
| | Cancelling a running week program Creating a new week program | |
| | 0.3.1 Section handling | |
| | 0.3.2 Temperature setpoint | |
| | 0.3.3 Day of the week | |
| - | 0.3.4 Time of the day | |
| | Activity of the shift-point Switching on or off the optional zero-voltage relay outputs | |
| | 0.3.7 Calling up the next parameter | |
| | 0.3.8 Saving the week program and leaving the program editor | |
| 10.4 | Deleting a week program | 77 |
| 11. | KEY LOCK | 78 |
| 11 1 | Directly activating the key lock function | 79 |
| 11.2 | Automatic key lock | |
| 11.3 | Changing the password for unlocking the key lock | |
| 12. | GENERAL CONTROLLER SETTINGS | |
| | | |
| 12.1 12.2 | Setup wizard Date and time settings | |
| 12.2 | Selecting the menu language of the T4.12 controller | |
| 12.4 | Setting display brightness | |
| 12.5 | Changing the temperature unit | 86 |
| 12.6 | Defining the data recording rate | |
| 12.7 12.8 | Factory reset | |
| 12.8 | Network configuration Display of the entire network configuration | |
| | Display and entry of the chamber configuration – for service purpose | |
| 13. | DATA TRANSFER VIA USB INTERFACE | |
| | | |
| 13.1 13.2 | Exporting data to USB drive Importing data from USB drive | |
| | | |

| 14. | NOTIFICATIONS AND ALARMS | 94 |
|--------------|--|-------|
| 14.1 | Notifications overview | - |
| | Alarms overview Alarm status | |
| | Confirming a "set" alarm. | |
| | Alarm configuration and overview | |
| | .5.2 History – list of all alarms | 98 |
| | .5.3 Activating, deactivating, and testing the alarm buzzer .5.4 Activating / deactivating all alarm functions | |
| 15. | EVENT LIST | |
| 16. | GRAPHICAL DISPLAY OF THE MEASURED VALUES | |
| 16.1 | Setting the sampling rate | |
| 16.2 | Defining the display range | |
| 16.3 | Selecting the parameters | 103 |
| 17. | TEMPERATURE SAFETY DEVICES | . 103 |
| 17.1 | Over temperature protective device (class 1) | 103 |
| | Overtemperature safety controller (temperature safety device class 3.1) | |
| 17 | .2.2 Setting the safety controller | 104 |
| | Over- and undertemperature safety controller (temperature safety device class 3.3) (option) 3.1 Safety controller modes | 108 |
| | .3.2 Setting the Over- and undertemperature safety controller | |
| 18. | DEFROSTING DURING REFRIGERATING OPERATION | 112 |
| 19. | OPTIONS | 113 |
| 19.1 | APT-COM™ 4 Multi Management Software (option) | 113 |
| 19.2 19.3 | Data logger kit (option) Zero-voltage relay alarm output (option) | |
| 19.3 19.4 | Water protected internal socket (option) | 113 |
| 19.5 19.6 | Object temperature display with flexible Pt 100 temperature sensor (option) Zero-voltage relay control outputs (may be available via BINDER INDIVIDUAL Customized | |
| | Solution) | |
| 19.7 | | |
| 20. | | |
| 20.1 20.2 | Cleaning Decontamination / chemical disinfection | |
| 21. | MAINTENANCE AND SERVICE, TROUBLESHOOTING, REPAIR, TESTING | .120 |
| 21.1 | General information, personnel qualification | 120 |
| 21.2 | Maintenance intervals, service | |
| 21.3 21.4 | Simple troubleshooting Sending the chamber back to BINDER GmbH | |
| 22. | DISPOSAL | |
| 22.1 | Disposal of the transport packing | 124 |
| 22.2 | Decommissioning | 125 |
| 22.3 22.4 | Disposal of the chamber in the Federal Republic of Germany Disposal of the chamber in the member states of the EU except for the Federal Republic of | 125 |
| | | 100 |
| 22.5 | Germany Disposal of the chamber in non-member states of the EU | |

| 23. | TECHNICAL DESCRIPTION | 128 |
|-------|---|-----|
| 23.1 | Factory calibration and adjustment | 128 |
| 23.2 | Over current protection | 128 |
| 23.3 | Definition of usable volume | 128 |
| | Technical data | |
| 23.5 | Equipment and Options (extract) | 131 |
| | Spare parts and accessories (extract) | |
| | KB / KB-UL 53 dimensions | |
| | KB / KB-UL 115 dimensions | |
| | KB / KB-UL 240 dimensions | |
| | KB / KB-UL 400 dimensions | |
| 23.11 | KB / KB-UL 720 dimensions | 137 |
| 24. | CERTIFICATES AND DECLARATIONS OF CONFORMITY | 138 |
| 24.1 | EU Declaration of Conformity | |
| | Certificate for the GS mark of conformity of the "Deutsche Gesetzliche Unfallversicherung | |
| | (German Social Accident Insurance) DGUV | 141 |
| 25. | PRODUCT REGISTRATION | 143 |
| 26. | CONTAMINATION CLEARANCE CERTIFICATE | 144 |
| 26.1 | For chambers located outside USA and Canada | 144 |
| | For chambers in USA and Canada | |

Dear customer,

For the correct operation of the chambers, it is important that you read this operating manual completely and carefully and observe all instructions as indicated. Observe the national regulations on minimum age of laboratory personnel. Failure to read, understand and follow the instructions may result in personal injury. It can also lead to damage to the chamber and/or poor equipment performance.

1. Safety

1.1 Personnel Qualification


The chamber must only be installed, tested, and started up by personnel qualified for assembly, startup, and operation of the chamber. Qualified personnel are persons whose professional education, knowledge, experience and knowledge of relevant standards allow them to assess, carry out, and identify any potential hazards in the work assigned to them. They must have been trained and instructed, and be authorized, to work on the chamber.

The chamber should only be operated by laboratory personnel especially trained for this purpose and familiar with all precautionary measures required for working in a laboratory. Observe the national regulations on minimum age of laboratory personnel.

1.2 Operating manual

This operating manual is part of the components of delivery. Always keep it handy for reference in the vicinity of the chamber. If selling the unit, hand over the operating manual to the purchaser.

To avoid injuries and damage observe the safety instructions of the operating manual. Failure to follow instructions and safety precautions can lead to significant risks.

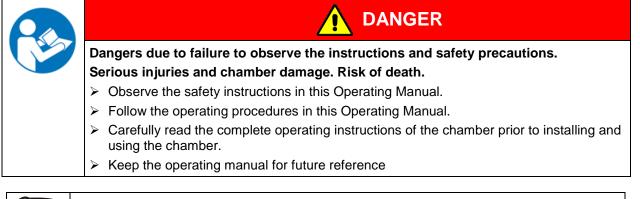

Make sure that all persons who use the chamber and its associated work equipment have read and understood the Operating Manual.

This Operating Manual is supplemented and updated as needed. Always use the most recent version of the Operating Manual. When in doubt, call the BINDER Service Hotline for information on the up-to-date-ness and validity of this Operating Manual.

1.3 Legal considerations

This operating manual is for informational purposes only. It contains information for correct and safe installing, start-up, operation, decommissioning, cleaning and maintenance of the product. Note: the contents and the product described are subject to change without notice.



Understanding and observing the instructions in this operating manual are prerequisites for hazard-free use and safety during operation and maintenance. Images are to provide basic understanding. They may deviate from the actual version of the chamber. The actual scope of delivery can, due to optional or special design, or due to recent technical changes, deviate from the information and illustrations in these instructions this operating manual. In no event shall BINDER be held liable for any damages, direct or incidental arising out of or related to the use of this manual.

This operating manual cannot cover all conceivable applications. If you would like additional information, or if special problems arise that are not sufficiently addressed in this manual, please ask your dealer or contact us directly, e.g. by phone at the number located on page one of this manual

Furthermore, we emphasize that the contents of this operating manual are not part of an earlier or existing agreement, description, or legal relationship, nor do they modify such a relationship. All obligations on the part of BINDER derive from the respective purchase contract, which also contains the entire and exclusively valid statement of warranty administration and the general terms and conditions, as well as the legal regulations valid at the time the contract is concluded. The statements in this manual neither augment nor restrict the contractual warranty provisions.

1.4 Structure of the safety instructions

In this operating manual, the following safety definitions and symbols indicate dangerous situations following the harmonization of ISO 3864-2 and ANSI Z535.6.

1.4.1 Signal word panel

Depending on the probability of serious consequences, potential dangers are identified with a signal word, the corresponding safety color, and if appropriate, the safety alert symbol.



WARNING

Indicates an imminently hazardous situation that, if not avoided, will result in death or serious (irreversible) injury.

Indicates a potentially hazardous situation which, if not avoided, could result in death or serious (irreversible) injury.



Indicates a potentially hazardous situation which, if not avoided, may result in moderate or minor (reversible) injury.

NOTICE

Indicates a potentially hazardous situation which, if not avoided, may result in damage to the product and/or its functions or of a property in its proximity.

1.4.2 Safety alert symbol



Use of the safety alert symbol indicates a risk of injury.

Observe all measures that are marked with the safety alert symbol in order to avoid death or injury.

1.4.3 Pictograms

| Warning signs | | | |
|---------------------------------|-------------------------|------------------------|---------------------------|
| Electrical hazard | Hot surface | Explosive atmosphere | Stability hazard |
| Lifting hazard | Risk of corrosion and / | Harmful substances | Biohazard |
| Pollution Hazard | or chemical burns | | |
| Mandatory action signs | 1 | 1 | |
| Mandatory regulation | Read operating | Disconnect the power | Lift with several persons |
| | instructions | plug | |
| | | | |
| Lift with mechanical assistance | Environment protection | Wear protective gloves | Wear safety goggles |
| | Environment protection | Wear protective gloves | Wear safety goggles |
| assistance | Environment protection | Wear protective gloves | Wear safety goggles |

1.4.4 Word message panel structure

Type / cause of hazard.

33

Possible consequences.

- $\ensuremath{\varnothing}$ Instruction how to avoid the hazard: prohibition
- > Instruction how to avoid the hazard: mandatory action

Observe all other notes and information not necessarily emphasized in the same way, in order to avoid disruptions that could result in direct or indirect injury or property damage.



1.5 Localization / position of safety labels on the chamber

The following labels are located on the chamber:

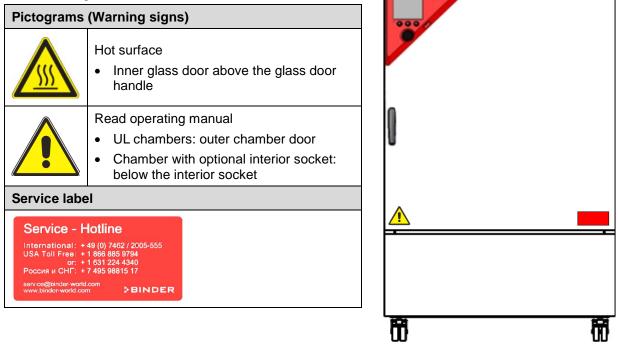


Figure 1: Position of labels on the chamber (example KB 240-UL)



Keep safety labels complete and legible.

Replace safety labels that are no longer legible. Contact BINDER Service for these replacements.

1.6 Type plate

The type plate sticks to the left side of the chamber, bottom right-hand.


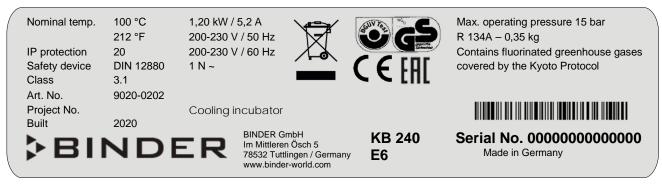


Figure 2: Type plate (example of KB 240 regular chamber)



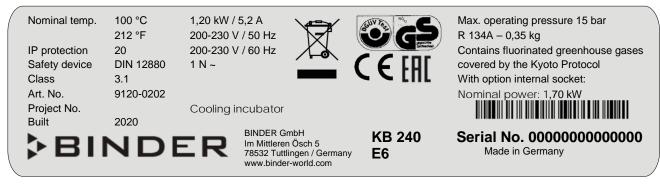

| Figure 3: Type | e plate (example of | f KB 240 optional cl | namber) |
|----------------|---------------------|----------------------|---------|
|----------------|---------------------|----------------------|---------|

| Indications of the | type plate (example) | Information | |
|---|---|---|--|
| BINDER | | Manufacturer: BINDER GmbH | |
| KB 240 | | Model designation | |
| Cooling incubator | | Device name | |
| Serial No. | Serial No. 000000000000 Serial no. of the chamber | | |
| Built | 2020 | Year of construction | |
| Nominal tempera- ture | 100 °C 212 °F | Nominal temperature | |
| IP protection | 20 | IP type of protection acc. to standard EN 60529 | |
| Temp. safety de- vice | DIN 12880 | Temperature safety device acc. to standard DIN 12880:2007 | |
| Class | 3.1 | Class of temperature safety device | |
| Art. No. | 9020-0202 | Art. no. of the chamber | |
| Project No. | | Optional: Special application acc. to project no. | |
| 1,20 kW | | Nominal power | |
| 5,2 A | | Nominal current | |
| 200-230 V / 50 Hz | | Nominal voltage range +/-10% | |
| 200-230 V / 60 Hz | | at the indicated power frequency | |
| 1 N ~ | | Current type | |
| Max. operating pre | ssure 15 bar | Max. operating pressure in the refrigerating system | |
| R 134A - 0,35 kg | | Refrigerant type and max. filling weight | |
| Contains fluorinated greenhouse gases covered by the Kyoto Protocol | | Contains fluorinated greenhouse gases covered by the Kyoto Protocol | |
| With option internal socket: Nominal power: 1,70 kW | | With option internal socket: increased total nominal power | |

| Symbol on the type plate | Information |
|--------------------------|--|
| CE | CE conformity marking |
| | Electrical and electronic equipment manufactured / placed on the market in the EU after 13 August 2005 and to be disposed of in separate collection according to Directive 2012/19/EU on waste electrical and electronic equipment (WEEE). |
| | GS mark of conformity of the "Deutsche Gesetzliche Unfallversi- cherung e.V. (DGUV), Prüf- und Zertifizierungsstelle Nahrungsmittel und Verpackung im DGUV Test" (German Social Accident Insur- ance (DGUV), Testing and Certification Body for Foodstuffs and Packaging Industry in DGUV Test). (Not valid for UL chambers) |




| Symbol on the type plate | Information | |
|--------------------------------|---|--|
| EHC | The chamber is certified according to Customs Union Technical Regulation (CU TR) for the Eurasian Economic Union (Russia, Belarus, Armenia, Kazakhstan Kyrgyzstan). | |
| | The chamber is certified by Underwriters Laboratories Inc. [®] according to the following standards: | |
| (KB-UL only) | • UL 61010-1, 3 rd Edition, 2012-05, Rev. 2015-07 | |
| LISTED LABORATORY EQUIPMENT | • CAN/CSA-C22.2 No. 61010-1, 3 rd Edition, 2012-05, Rev. 2015-07 | |
| 43KM | • IEC/EN 61010-1:2014, 3 rd Edition | |

1.7 General safety instructions on installing and operating the chambers


With regard to operating the chambers and to the installation location, please observe the local and national regulations relevant for your country (for Germany: DGUV guidelines 213-850 on safe working in laboratories, issued by the employers' liability insurance association).

BINDER GmbH is only responsible for the safety features of the chamber provided skilled electricians or qualified personnel authorized by BINDER perform all maintenance and repair, and if components relating to chamber safety are replaced in the event of failure with original spare parts.


To operate the chamber, use only original BINDER accessories or accessories from third-party suppliers authorized by BINDER. The user is responsible for any risk caused by using unauthorized accessories.

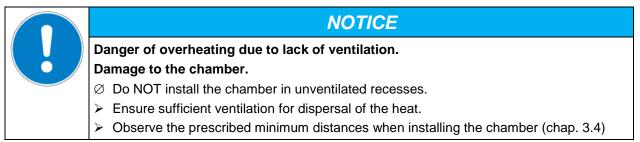

Do not install or operate the chambers in hazardous locations.

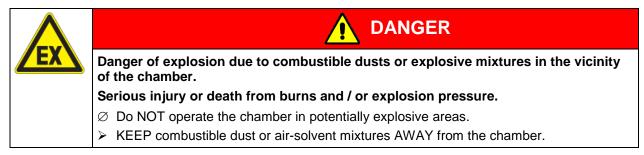


The chambers do not dispose of any measures of explosion protection.

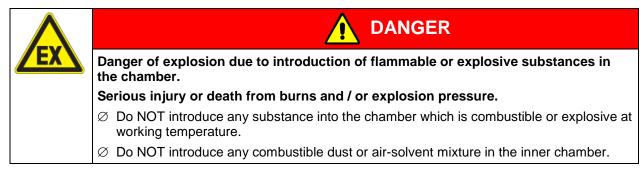




Any solvent contained in the charging material must not be explosive or inflammable. I.e., irrespective of the solvent concentration in the steam room, NO explosive mixture with air must form. The temperature inside the chamber must lie below the flash point or below the sublimation point of the charging material. Familiarize yourself with the physical and chemical properties of the charging material, as well as the contained moisture constituent and its behavior with the addition of heat energy.

Familiarize yourself with any potential health risks caused by the charging material, the contained moisture constituent or by reaction products which may arise during the temperature process. Take adequate measures to exclude such risks prior to putting the cooling incubator into operation.

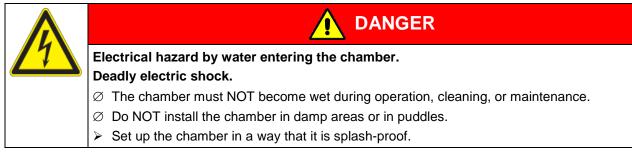

The chambers were produced in accordance with VDE regulations and were routinely tested in accordance to VDE 0411-1 (IEC 61010-1).

During and shortly after operation, the temperature of the inner surfaces almost equals the set-point. The glass doors, the glass door handles, and the inner chamber will become hot during operation.

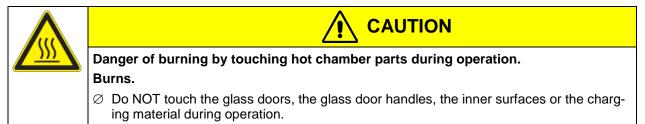





Danger of injury and damages by the chamber tipping over or breakaway of the protruding lower housing cover.

Injuries and damage to the chamber and the loading material

arnothing Do NOT load the lower housing cover with heavy objects while the chamber door is open and do NOT climb on it.

1.8 Intended use

Following the instructions in this operating manual and conducting regular maintenance work (chap. 21.2) are part of the intended use.

Any use of the chambers that does not comply with the requirements specified in this Operating Manual shall be considered improper use.

Other applications than those described in this chapter are not approved.

Use


Cooling incubators KB / KB-UL are suitable for exact conditioning of harmless materials. Because of their precise temperature accuracy these devices are especially useful for cultivation of microorganisms with a narrow temperature optimum in a range of 4 °C / 39.2 °F to 37 °C / 98.6 °F. Main fields of application are tests of long-term storage (e.g. at 4 °C / 39.2 °F), refrigerated incubation between 20 °C / 68 °F and 25 °C / 77 °F and incubation at 37 °C / 98.6 °F (also with additional introduction of heat) or with alternating temperatures (e.g. 37 °C / 98.6 °F and 4 °C / 39.2 °F).

Requirements for the chamber load


Any solvent must not be explosive and flammable. A mixture of any component of the charging material with air must NOT be explosive. The operating temperature must lie below the flash point or below the sublimation point of the charging material. Any component of the charging material must NOT be able to release toxic gases.

The charging material shall not contain any corrosive ingredients that may damage the machine components made of stainless steel, aluminum, and copper. Such ingredients include in particular acids and halides. Any corrosive damage caused by such ingredients is excluded from liability by BINDER GmbH.

The chamber does not dispose of any measures of explosion protection.

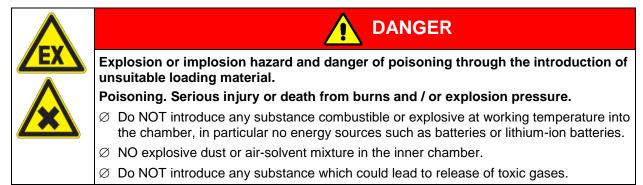

Contamination of the chamber by toxic, infectious or radioactive substances must be prevented

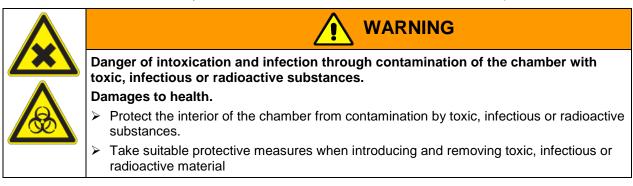

In case of foreseeable use of the chamber there is no risk for the user through the integration of the chamber into systems or by special environmental or operating conditions in the sense of EN 61010-1:2010. For this, the intended use of the chamber and all its connections must be observed.

Medical devices

The chambers are not classified as medical devices as defined by the Medical Device Directive 93/42/EEC.



Due to the special demands of the Medical Device Directive (MDD), these chambers are not qualified for sterilization of medical devices as defined by the directive 93/42/EWG.

Personnel Requirements

Only trained personnel with knowledge of the Operating Manual can set up and install the chamber, start it up, operate, clean, and take it out of operation. Service and repairs call for further technical requirements (e.g. electrical know-how), as well as knowledge of the service manual.

Installation site requirements

The chambers are designed for setting up inside a building (indoor use).

The requirements described in the Operating Manual for installation site and ambient conditions (Chap. 3.4) must be met.

WARNING: If customer should use a BINDER chamber running in non-supervised continuous operation, we strongly recommend in case of inclusion of irrecoverable specimen or samples to split such specimen or samples and store them in at least two chambers, if this is feasible.

1.9 Foreseeable Misuse

Other applications than those described in chap. 1.8 are not approved.

This expressly includes the following misuses (the list is not exhaustive), which pose risks despite the inherently safe construction and existing technical safety equipment:

- Non-observance of Operating Manual
- Non-observance of information and warnings on the chamber (e.g. control unit messages, safety identifiers, warning signals)
- Installation, startup, operation, maintenance and repair by untrained, insufficiently qualified, or unauthorized personnel
- Missed or delayed maintenance and testing
- Non-observance of traces of wear and tear
- Insertion of materials excluded or not permitted by this Operating Manual.
- Non-compliance with the admissible parameters for processing the respective material.
- Installation, testing, service or repair in the presence of solvents
- Installation of replacement parts and use of accessories and operating resources not specified and authorized by the manufacturer
- Installation, startup, operation, maintenance or repair of the chamber in absence of operating instructions
- Bypassing or changing protective systems, operation of the chamber without the designated protective systems
- Non-observance of messages regarding cleaning and disinfection of the chamber.
- Spilling water or cleaning agent on the chamber, water penetrating into the chamber during operation, cleaning or maintenance.

- Cleaning activity while the chamber is turned on.
- Operation of the chamber with a damaged housing or damaged power cord
- Continued operation of the chamber during an obvious malfunction
- Insertion of objects, particularly metallic objects, in louvers or other openings or slots on the chamber
- Human error (e.g. insufficient experience, qualification, stress, exhaustion, laziness)

To prevent these and other risks from incorrect operation, the operator shall issue operating instructions. Standard operating procedures (SOPs) are recommended.

1.10 Residual Risks

The unavoidable design features of a chamber, as well as its proper field of application, can also pose risks, even during correct operation. These residual risks include hazards which, despite the inherently safe design, existing technical protective equipment, safety precautions and supplementary protective measures, cannot be ruled out.

Messages on the chamber and in the Operating Manual warn of residual risks. The consequences of these residual risks and the measures required to prevent them are listed in the Operating Manual. Moreover, the operator must take measures to minimize hazards from unavoidable residual risks. This includes, in particular, issuing operating instructions.

The following list summarizes the hazards against which this Operating Manual and the Service Manual warn, and specifies protective measures at the appropriate spots:

Unpacking, Transport, Installation

- Sliding or tilting the chamber
- Setup of the chamber in unauthorized areas
- · Installation of a damaged chamber
- Installation of a chamber with damaged power cord
- Inappropriate site of installation
- Missing protective conductor connection

Normal operation

- Assembly errors
- · Contact with hot surfaces on the housing
- Contact with hot surfaces in the interior and inside of doors
- Emission of non-ionizing radiation from electrical operating resources
- Contact with live parts in normal state

Cleaning and Decontamination

- Penetration of water into the chamber
- Inappropriate cleaning and decontamination agents
- Enclosure of persons in the interior

Malfunction and Damage

- Continued operation of the chamber during an obvious malfunction or outage of the heating or cooling system
- Contact with live parts during error status
- Operation of a unit with damaged power cord

Maintenance

- Maintenance work on live parts.
- Execution of maintenance work by untrained/insufficiently qualified personnel
- · Electrical safety analysis during annual maintenance not performed

Trouble-shooting and Repairs

- Non-observance of warning messages in the Service Manual
- Trouble-shooting of live parts without specified safety measures
- Absence of a plausibility check to rule out erroneous inscription of electrical components
- Performance of repair work by untrained/insufficiently qualified personnel
- Inappropriate repairs which do not meet the quality standard specified by BINDER
- Use of replacement parts other than BINDER original replacement parts
- Electrical safety analysis not performed after repairs

1.11 Operating instructions

Depending on the application and location of the chamber, the operator of the chamber must provide the relevant information for safe operation of the chamber in a set of operating instructions.

(F)

Keep these operating instructions with the chamber at all times in a place where they are clearly visible. They must be comprehensible and written in the language of the employees.

1.12 Measures to prevent accidents

The operator of the chamber must observe the following rule: "Betreiben von Arbeitsmitteln. Betreiben von Kälteanlagen, Wärmepumpen und Kühleinrichtungen" (Operation of work equipment. Operation of refrigeration systems, heat pumps and refrigeration equipment) (GUV-R 500 chap. 2.35) (for Germany).

The manufacturer took the following measures to prevent ignition and explosions:

• Indications on the type plate

See operating manual chap. 1.6.

• Operating manual

An operating manual is available for each chamber.

• Overtemperature monitoring

The chamber is equipped with a temperature display, which can be read from outside.

The chamber is equipped with an additional safety controller (temperature safety device class 3.1 acc. to DIN 12880:2007). Visual and audible (buzzer) signals indicate temperature exceeding.

• Safety, measurement, and control equipment

The safety, measuring, and control equipment is easily accessible.

• Electrostatic charge

The interior parts are grounded.

• Non-ionizing radiation

Non-ionizing radiation is not intentionally produced, but released only for technical reasons by electrical equipment (e.g. electric motors, power cables, solenoids). The machine has no permanent magnets. If persons with active implants (e.g. pacemakers, defibrillators) keep a safe distance (distance of field source to implant) of 30 cm, an influence of these implants can be excluded with high probability.

• Protection against touchable surfaces

Tested according to EN ISO 13732-1:2008.

Floors

See operating manual chap. 3.4 for correct installation

• Cleaning

See operating manual chap. 20.

• Examinations

The chamber has been inspected by the "Deutsche Gesetzliche Unfallversicherung e.V. (DGUV) (German Social Accident Insurance (DGUV)" (German Social Accident Insurance (DGUV), Testing and Certification Body for Foodstuffs and Packaging Industry in DGUV Test) and bears the GS mark. (Not valid for UL chambers)

UL units only: The chamber is certified by Underwriters Laboratories Inc.[®] according to the following standards: UL 61010-1, 3rd Edition, 2012-05, Rev. 2015-07; CAN/CSA-C22.2 No. 61010-1, 3rd Edition, 2012-05, Rev. 2015-07; IEC/EN 61010-1:2014, 3rd Edition

2. Chamber description

A high level of precision, reliability, and safety for all growth parameters ensures optimum incubation conditions. Moreover, the KB / KB-UL cooling incubator is designed for maximum usability – even in continuous operation year after year. It fulfills all technical and application-specific requirements arising in experimentation such as in the areas of biotechnology, medicine, the nutrition industry, pharmaceutical and cosmetics industries, botany, and zoology.

Two important temperature technologies have been combined to achieve perfect temperature control. The specially developed DCT[™] refrigerating system, a direct refrigerating process, in conjunction with the APT.line[™] preheating chamber technology, satisfies the unique prerequisites for attaining highly-precise temperature control and particularly short recovery times after opening the door.

The refrigerating system is distinguished by direct, precise, and rapid temperature conduction. Large-area labyrinth evaporator plates directly conduct the cold to the atmosphere of the working space.

The APT.line[™] preheating chamber system ensures high level of spatial and time-based temperature precision, thanks to the direct and distributed air circulation into the interior. This is especially important for maintaining temperatures – especially with full chambers – and for rapid restoration of optimum growth conditions after opening the door. The inner glass door ensures that the temperature remains constant when observing the incubation process. The fan supports exact attainment and maintenance of the desired temperature accuracy. The fan speed is digitally adjustable. The heating and refrigerating systems are microprocessor regulated to a tenth of a degree. In addition, the chamber provides almost unlimited possibilities for adaptation to individual customer requirements based upon extensive programming options and on the week program timer and real time clock of the controller.

The KB / KB-UL 240 / 400 / 720 (E6) chambers are equipped with a door heating.

All chamber functions are easy and comfortable to use thanks to their clear arrangement. Major features are easy cleaning of all chamber parts and avoidance of undesired contamination.


Material: The inner chamber, the pre-heating chamber and the interior side of the doors are all made of stainless steel V2A (German material no. 1.4301, US equivalent AISI 304). The housing is RAL 7035 pow-der-coated. All corners and edges are also completely coated.

The chamber comes equipped with an Ethernet interface for computer communication, e.g. via the APT-COM[™] 4 Multi Management Software (option, chap. 19.1). For further options, see chap. 23.5.

The KB / KB-UL 240 / 400 / 720 (E6) chambers are equipped with four castors. Both front castors can easily be locked via the attached brakes.

Temperature range at ambient temperature of +22 °C +/- 3°C / 71.6 °F +/- 5.4 °F: -10 °C / 14 °F up to +100 °C / 212 °F

2.1 Chamber overview

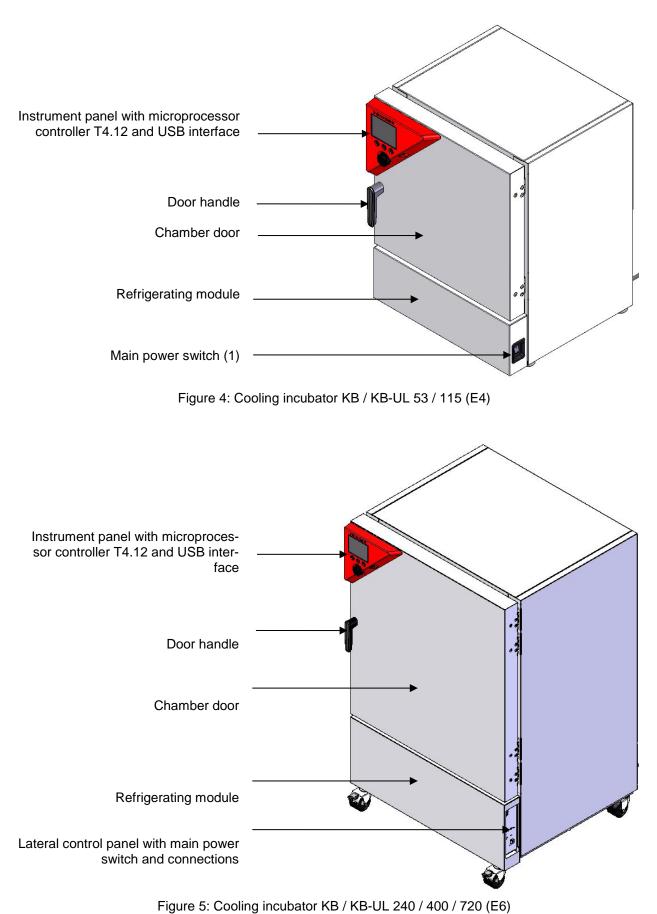



2.2 Instrument panel

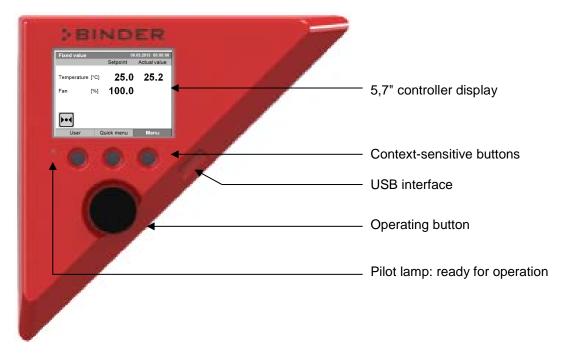


Figure 6: Instrument panel with microprocessor controller T4.12 and USB interface

2.3 Lateral control panel with main power switch and optional equipment – KB / KB-UL 240 / 400 / 720 (E6)

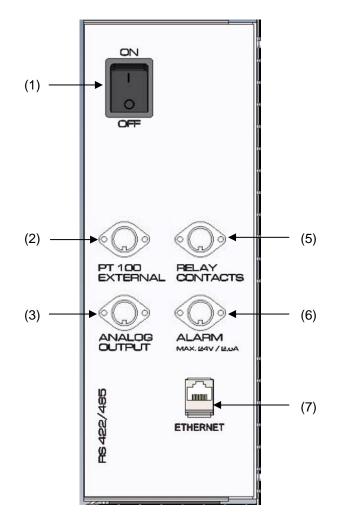


Figure 7: Lateral control panel at the right side of the refrigerating machine, with options

- (1) Main power switch
- (2) DIN-socket for additional Pt 100 temperature sensor (may be available via BINDER INDIVIDUAL Customized Solutions)
- (3) DIN socket for analog output 4-20 mA (option)
- (4) (not used)
- (5) DIN-socket for zero-voltage relay control outputs (may be available via BINDER INDIVIDUAL Customized Solutions)
- (6) DIN-socket for zero-voltage relay alarm output (option)
- (7) Ethernet interface for computer communication

2.4 Chamber rear – KB / KB-UL 53 / 115 (E4)

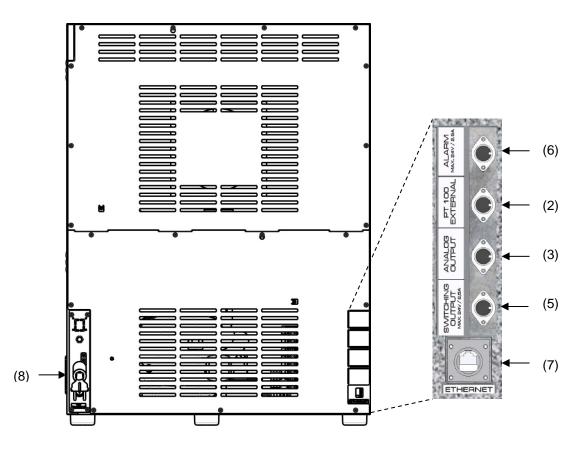


Figure 8: Chamber rear with position of options (example)

- (2) DIN-socket for additional Pt 100 temperature sensor (may be available via BINDER INDIVIDUAL Customized Solutions)
- (3) DIN socket for analog output 4-20 mA (option)
- (4) (not used)
- (5) DIN-socket for zero-voltage relay control outputs (may be available via BINDER INDIVIDUAL Customized Solutions)
- (6) DIN-socket for zero-voltage relay alarm output (option)
- (7) Ethernet interface for computer communication
- (8) Power cable


3. Completeness of delivery, transportation, storage, and installation

3.1 Unpacking, and checking equipment and completeness of delivery

After unpacking, please check the chamber and its optional accessories, if any, based on the delivery receipt for completeness and for transportation damage. Inform the carrier immediately if transportation damage has occurred.

The final tests of the manufacturer may have caused traces of the shelves on the inner surfaces. This has no impact on the function and performance of the chamber.

Please remove any transportation protection devices and adhesives in/on the chamber and on the doors and remove the operating manuals and accessory equipment.

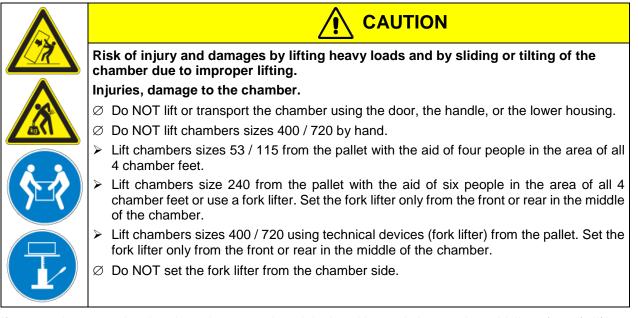

If you need to return the chamber, please use the original packing and observe the guidelines for safe lifting and transportation (chap. 3.2).

For disposal of the transport packing, see chap. 22.1.


Note on second-hand chambers (Ex-Demo-Units):

Second-hand chambers are chambers that were used for a short time for tests or exhibitions. They are thoroughly tested before resale. BINDER ensures that the chamber is technically sound and will work flaw-lessly.

Second-hand chambers are marked with a sticker on the chamber door. Please remove the sticker before commissioning the chamber.

3.2 Guidelines for safe lifting and transportation

The front castors of the chambers sizes 240 / 400 / 720 can be blocked by brakes. Please move the chambers with castors only when empty and on an even surface, otherwise the castors may be damaged. After operation, please observe the guidelines for temporarily decommissioning the chamber (chap. 22.2).

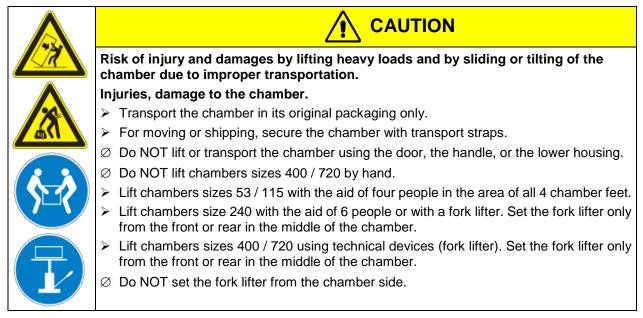

• Permissible ambient temperature range during transport: -10 °C / 14°F to +60 °C / 140°F.

You can order transport packing for moving or shipping purposes from BINDER service.

3.3 Storage

Intermediate storage of the chamber is possible in a closed and dry room. Observe the guidelines for temporary decommissioning (chap. 22.2).

- Permissible ambient temperature range during storage: -10 °C / 14°F to +60 °C / 140°F.
- Permissible ambient humidity: max. 70 % r.H., non-condensing

When after storage in a cold location you transfer the chamber to its warmer installation site, condensation may form. Before start-up, wait at least one hour until the chamber has attained ambient temperature and is completely dry.

3.4 Location of installation and ambient conditions

Set up the chamber on a flat, even surface, free from vibration, in a well-ventilated, dry location and align it using a spirit level. The site of installation must be capable of supporting the chamber's weight (see technical data, chap. 23.4). The chambers are designed for setting up inside a building (indoor use).

| NOTICE |
|---|
| Danger of overheating due to lack of ventilation. Damage to the chamber. |
| \varnothing Do NOT install the chamber in unventilated recesses. |
| Ensure sufficient ventilation for dispersal of the heat. |
| Observe the prescribed minimum distances when installing the chamber. |

Do not install or operate the chamber in potentially explosive areas.



Ambient conditions

• Permissible ambient temperature range during operation: +18 °C / 64.4 °F to +32 °C / 89.6 °F. At elevated ambient temperature values, fluctuations in temperature can occur.



The ambient temperature should not be substantially higher than the indicated ambient temperature of +22 °C +/- 3 °C / 71.6 °F +/- 5.4 °F to which the specified technical data relates. Deviations from the indicated data are possible for other ambient conditions.



With each degree of ambient temperature > +25 °C / 77°F, the refrigeration power decreases by 1.5 K.

• Permissible ambient humidity: 70 % r.H. max., non-condensing.

When operating the chamber at temperature set-points below ambient temperature, high ambient humidity may lead to condensation on the chamber.

• Installation height: max. 2000 m / 6562 ft. above sea level.

Minimum distances

- When placing several chambers of the same size side by side, maintain a minimum distance of 250 mm / 9.84 *in* between each chamber.
- Wall distances: rear 100 mm / 3.9 in, sides 160 mm / 6.29 in.
- Spacing above the chamber of at least 100 mm / 3.9 in must also be accounted for.

Stacking

Two KB / KB-UL 53 / 115 (E4) chambers of the same size can be piled on top of each other. For this purpose place rubber pads under every foot of the upper chamber to prevent the device from slipping.



KB / KB-UL 240 / 400 / 720 (E6) chambers must NOT be stacked.



Other requirements

To completely separate the chamber from the power supply, you must disconnect the power plug. Install the chamber in a way that the power plug is easily accessible and can be easily pulled in case of danger.

For the user there is no risk of temporary overvoltages in the sense of EN 61010-1:2010.

With an increased amount of dust in the ambient air, clean the condenser fan (by suction or blowing) several times a year.

Avoid any conductive dust in the ambiance according to the chamber layout complying with pollution degree 2 (IEC 61010-1).

4. Installation of the chamber

4.1 Spacers for wall distance – KB / KB-UL 240 / 400 / 720 (E6)

Please fix both spacers with the supplied screws at the chamber rear. This serves to ensure the prescribed minimum distance to the wall of 100 mm / 3.94 in.



Figure 9: Spacer for rear wall distance

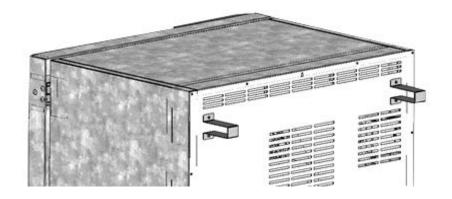


Figure 10: Rear KB / KB-UL 240 / 400 / 720 (E6) with mounted spacers

4.2 Mounting the flexible tilt protection kit

Use the supplied flexible tilt protection kit in addition to the spacers for wall distance (chap. 4.1). This will prevent the chamber from tilting when the door is open.



NOTICE

Danger of damages caused by tilting of the chamber when the door is open. Damage to the chamber.

Fix the chamber to a wall with supplied flexible tilt protection kit

Scope of delivery:

• 4 Torx screws (spare parts)

 \geq

- 4 tilt protection holders
- 4 securing straps (2 spare parts)

Mounting on device side:

- Remove two screws on the upper part of the rear wall (a)
- Fix two of the supplied tilt protection holders, each centrally with one of these screws (b).

Mounting on wall side

• Fix two of the supplied tilt protection holders in the appropriate distance, each with two screws Ø 6mm suitable for the wall (c)

Connection with the securing straps

• On each side, thread one of the supplied securing straps through the provided slots of a device side and a wall side tilt protection holder

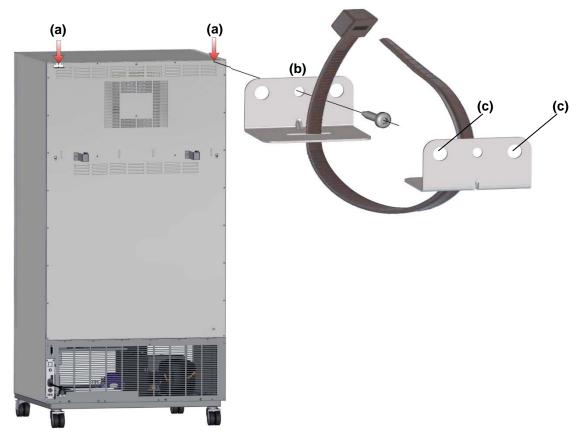


Figure 11: Rear chamber (example KB 400) and mounting the flexible tilt protection kit

4.3 Electrical connection

The chambers are supplied ready for connection. They come with a fixed power connection cable that has a length of at least 1800 mm / 5.9 ft and are protected with a chamber-protection against over current, the chambers KB 53-UL / KB 115-UL in addition with an internal fuse.

| Model | Power plug | Nominal voltage +/- 10% at the indicated power frequency | Current type | Chamber fuse |
|---|------------------|---|-----------------|-----------------|
| KB 53 (E4) KB 115 (E4) | Shock-proof plug | 230 V at 50 Hz | 1N~ | 10 A |
| KB 240 (E6) KB 400 (E6) KB 720 (E6) | Shock-proof plug | 200-230 V at 50 Hz 200-230 V at 60 Hz | 1N~ | 16 A |
| KB 53-UL (E4) KB 115-UL (E4) | NEMA 5-15 | 115 V at 60 Hz | 1N~ | 12,5 A |
| KB 240-UL (E6) KB 400-UL (E6) | NEMA 5-20P | 100-120 V at 50 Hz 100-120 V at 60 Hz | 1N~ | 16 A |
| KB 720-UL (E6) | NEMA 6-20P | 200-240 V at 50 Hz 200-240 V at 60 Hz | 2 ~ | 16 A |

• The domestic socket must also provide a protective conductor. Make sure that the connection of the protective conductor of the domestic installations to the chamber's protective conductor meets the latest technology. The protective conductors of the socket and plug must be compatible!



Electrical hazard due to missing protective conductor connection.

- Deadly electric shock.
- Make sure that the chamber's power plug and the power socket match and securely connect the electrical protective conductors of the chamber and the house installation.
- Only use original connection cables from BINDER according to the above specification.
- Prior to connection and start-up, check the power supply voltage. Compare the values to the specified data located on the chamber's type plate (left chamber side, bottom right-hand, see chap. 1.6).



NOTICE

Danger of incorrect power supply voltage due to improper connection. Damage to the chamber.

- Check the power supply voltage before connection and start-up.
- > Compare the power supply voltage with the data indicated on the type plate.
- When connecting, please observe the regulations specified by the local electricity supply company as well as the local or national electrical regulations (VDE directives for Germany).
- Observe a sufficient current protection according to the number of devices that you want to operate. We recommend the use of a residual current circuit breaker.
- Pollution degree (acc. to IEC 61010-1): 2
- Over-voltage category (acc. to IEC 61010-1): II

See also electrical data (chap. 23.4).





To completely separate the chamber from the power supply, you must disconnect the power plug. Install the chamber in a way that the power plug is easily accessible and can be easily pulled in case of danger.

5. Start up

After connecting the electrical supply (chap. 4.1), turn on the chamber by the main power switch (1). The pilot lamp shows the chamber is ready for operation.

Observe a delay time of about 30s between turning Off and On again. Otherwise an initialization problem may occur.

Note that the chamber is in stand-by mode when the main power switch has been turned on and yet the controller display is dark. Turn on the chamber by pressing any controller button.

Warming chambers may release odors in the first few days after commissioning. This is not a quality defect. To reduce odors quickly we recommend heating up the chamber to its nominal temperature for one day and in a well-ventilated location.

WARNING: If customer should use a BINDER chamber running in non-supervised continuous operation, we strongly recommend in case of inclusion of irrecoverable specimen or samples to split such specimen or samples and store them in at least two chambers, if this is feasible.

6. Functional overview of the T4.12 chamber controller

The T4.12 chamber controller controls the temperature (range: -10 °C / 14 °F up to +100 °C / 212 °F) and the fan speed (range: 40 % up to 100 %) inside the chamber. You can enter the desired set point values in fixed value operating mode or in program mode in the display controller. The controller also offers a week program function and various notifications and alarm messages with visual and audible indication, a trace file and remote alarms via e-mail. You can enter values or programs directly at the controller keypad or using the APT-COMTM 4 Multi Management Software (option, chap. 19.1) specially developed by BINDER.

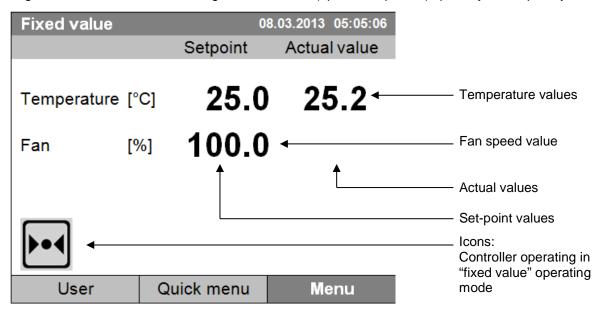


Figure 12: T4.12 microprocessor controller, initial view in "fixed value" controller mode (sample values)

6.1 Menu structure

| Fixed value | ÷ | 08. | .03.2015 05:05:06 |
|-------------|--------|-----------|-------------------|
| | | Setpoint | Actual value |
| Temperature | e [°C] | 25.0 | 25.2 |
| Fan | [%] | 100.0 | |
| | | | |
| | | | |
| ▶•◀ | | | |
| User | C | uick menu | Menu |

From the Initial view you have access to different menus using the menu buttons "User", "Quick menu", or "Menu". From there you can access the desired control functions. To do this, select the function by turning the operating button and press the operating button to confirm the selection.

In any menu, you can return to the previous display pressing the "Close" button or to the initial view with the "Home" button.

Depending on the logged-in user or administrator, the available menu functions may vary. These instructions present the functions which are available to the logged-in administrator.

6.1.1 General menu

The general menu provides access to all setting functions of the controller, a graphical display of the measured values, and the possibility to read and give out data via the USB interface. In addition, supporting functions like a settings wizard or a contact page are available.

| Fixed value | 08.03.2015 | 05:05:06 | |
|-------------------|------------|---------------|--------------|
| \ Menu | | | |
| Controller mode | | | |
| Event list | | | |
| Alarms | | | |
| Setpoints | | | |
| Safety controller | | | General menu |
| Programs | | | |
| Import/Export | | | |
| Settings | | $\overline{}$ | |
| | | · | |
| Close | Ho | ome | |

Turn the operating button to see additional menu items.

| Fixed value | 08.03.2015 05 | :05:06 | |
|------------------|---------------|--------|--|
| \ Menu | | | |
| Measurement cl | hart | | |
| Optional equipme | ent | | |
| Sensor adjustme | nt | | |
| Service contact | | | General menu (next page) |
| System informati | on | | ("Optional equipment" menu item is visible only with optional chamber equipment) |
| | | | |
| | | | |
| | | | |
| | _ | | |
| Close | | Home | |

| Controller mode | Switching between the operating modes "control off" or "fixed value", chap. 6.2.1 |
|--------------------|--|
| Event list | Display of status information and errors, chap. 15 |
| Alarms | Alarm settings, chap. 14.5 |
| Setpoints | Setpoint entry in "Fixed value" operating mode, chap. 8 |
| Safety controller | Setting the safety controller, chap. 17.2 and 17.3 |
| Programs | Time and week programs, chap. 9 and 10 |
| Import/Export | Data transfer via USB interface, chap. 13 |
| Settings | General controller settings, chap. 12 |
| Measurement chart | Graphical display of the measured values, chap. 16 |
| Optional equipment | Setting for optional equipment like interior socket, zero-voltage relay control outputs, alarm output, object temperature display, chap. 7 (menu item is visible only with optional chamber equipment) |
| Sensor adjustment | Adjustment menu for single-point and two-point adjustments (for Service purpose) |
| Service contact | Service information |
| System information | Chamber information (model, name, serial no., firmware etc.) |

6.1.2 Quick menu

| Fixed value | 08.03.2015 | 05:05:06 | |
|----------------------------|------------|----------|-----------------|
| \ Quick menu | | | |
| Measurement chart | | | |
| Active alarms | | | |
| Temperature setpoint | | | |
| Fan speed setpoint | | | "Ouisland and " |
| Safety controller setpoint | | | "Quick menu" |
| Time program | | | |
| Week program | | | |
| | | | |
| | | | |
| Close | Ho | me | |

The Quick menu provides fast access to frequently used functions.

| Measurement chart | Graphical display of the measured values, chap. 16 |
|----------------------------|---|
| Active alarms | Alarm settings, chap. 14.5 |
| Temperature setpoint | Temperature setpoint entry in "Fixed value" operating mode, chap. 8 |
| Fan speed setpoint | Fan speed setpoint entry in "Fixed value" operating mode, chap. 8 |
| Safety controller setpoint | Setting the safety controller setpoint, chap. 17.2 and 17.3 |
| Time program | Starting and cancelling a time program, chap. 9.1, 9.2 |
| Week program | Starting and cancelling a week program, chap. 10.1, 10.2 |

6.1.3 "User" menu

The user menu includes the key lock function and provides quick access to the event list. The key lock function serves to block the access to the controller. An overview of logon, logoff, and other events is given in the event list.

| Fixed value | 0 | 8.03.2015 | 05:05:06 | |
|-----------------|---|-----------|----------|-------------|
| \ User | | | | |
| Key lock | | | | |
| Show event list | | | | |
| | | | | |
| | | | | "User" menu |
| | | | | |
| | | | | |
| | | | | |
| | | | | |
| | l | | | |
| Close | | Но | me | |

| Key lock | Configuring the key lock function, chap. 11 |
|-----------------|---|
| Show event list | Displaying the event list, chap. 15 |

6.2 Operating modes

In the "**control off**" mode (chap. 6.2.1), the controller is non-functional and displays only the actual values. There is no heating or refrigeration. The temperature approximates the ambient value, the fan turns with 40 % speed.

You can enter the desired set point values in "**fixed value**" mode (chap. 8). The controller then operates as a fixed-point control, i.e., it reaches and maintains the defined temperature set-point until the next manual change.

The T4.12 program controller also permits running a **time program** (chap. 9) or a **week program** (chap. 10). You can program temperature cycles and define also the fan speed for each program section. The controller offers 52 time program places with up to 100 sections each. The week program mode offers 8 week program places with up to 30 shift points for each week program.

6.2.1 Activating the "control off" mode or change to "fixed value" operating mode

To select the "control off" or "fixed value" operating mode, go to Menu > Controller mode

| Fixed value | 0 | 8.03.2015 | 05:05:06 | |
|-------------------------------|---|-----------|----------|--|
| \ Menu | | | | |
| Controller mod | e | | | |
| Event list | | | | |
| Alarms | | | | |
| Setpoints | | | | General menu. |
| Safety controller | r | | | Select "Controller Mode" and press the operating button. |
| Programs | | | | and press the operating button. |
| Import/Export | | | | |
| Settings | | | | |
| | | | | |
| Close | | Ho | me | |
| | | | | |
| | | | | |
| Fixed value | | 8.03.2015 | 05:05:06 | |
| \ Controller m | | 8.03.2015 | 05:05:06 | |
| \ Controller m Fixed value | | 8.03.2015 | 05:05:06 | |
| \ Controller m | | 8.03.2015 | 05:05:06 | |
| \ Controller m Fixed value | | 8.03.2015 | 05:05:06 | Submenu "Controller Mode". |
| \ Controller m Fixed value | | 8.03.2015 | 05:05:06 | Select the desired controller mode |
| \ Controller m Fixed value | | 8.03.2015 | 05:05:06 | Select the desired controller mode "Control off" or "Fixed value" |
| \ Controller m Fixed value | | 8.03.2015 | 05:05:06 | Select the desired controller mode |
| \ Controller m Fixed value | | 8.03.2015 | 05:05:06 | Select the desired controller mode "Control off" or "Fixed value" |
| \ Controller m Fixed value | | 8.03.2015 | 05:05:06 | Select the desired controller mode "Control off" or "Fixed value" |
| \ Controller m Fixed value | | 8.03.2015 | 05:05:06 | Select the desired controller mode "Control off" or "Fixed value" |



| Control off \ Menu Controller mode Event list Alarms Setpoints Safety controller Programs Import/Export Settings | 08.03.2015 05:05:06 | General menu with controller mode "Control off". The controller mode "Fixed value" or "Control off" is indicated in the display headline. |
|---|-------------------------------------|---|
| Close Go back to the initial view with "H | Home łome". | |
| Control off | 08.03.2015 05:05:06 Actual value | |
| Temperature [°C] Fan [%] 40. | 25.2 0 | Initial view in "Control off" mode (sample picture). |
| User Quick menu | Menu | |

The controller is non-functional, i.e., there is no heating or refrigerating. The fan turns at 40 % speed.

| } | In the "Control off" mode, no program can be started. |
|----------|---|
| | |

6.3 Performance during and after power failure

During a power failure, all controller functions are shut down. The optional zero-voltage relay alarm output (chap. 19.3) is switched to alarm position for the whole duration of the power failure.

After the power returns, all functions return to the same status the chamber had before power failure. The controller continues to function in the original operating mode it was in previously before the power failure occurred.

• Performance after power failure in "fixed value" operation mode

All functions return to the same status the chamber had before power failure. The set-points are immediately resumed.

• Performance after power failure during time program operation

The program is resumed at the point where the interruption occurred with the latest set-points reached during the program run.

• Performance after power failure during week program operation

The week program continues with the values corresponding to the current time.

In the "Control off" controller mode, no program can be started.

If the temperature has dropped below the alarm limit during power failure, confirm the alarm with the RESET button as soon as the correct values are reached again (chap. 14.4).


In the Event List (chap. 15) the message "Power supply was interrupted." is shown.

6.4 Information

You access chamber information like the chamber type, serial no., firmware version etc. To display the system information, go to *Menu* > *System information*

| Fixed value | 08.03.2015 05 | 5:05:06 | |
|---------------------|-------------------|---------|---|
| \ System infor | mation | | |
| Chamber type: | KB | | |
| Chamber name: | KB_E6 | | |
| Serial number: (| 00-00000 | | |
| Special application | on number: 00-00 | 000 | Submenu "System information" (sample values). |
| Parameter versi | on: 511B-000A-0 | 005 | |
| Firmware versio | n (1): 521C-0001 | -0030 | |
| Firmware versio | n (2): 521B-0005- | -0028 | |
| | | | |
| | | | |
| Close | | Home | |

To display the BINDER Service contact data, go to *Menu* > Service contact

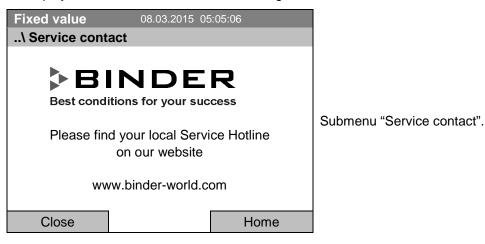


Further information windows are accessible under *Menu* > *Settings* > *Network settings* > *Show network settings* (chap. 12.9) and – for service purpose – under *Menu* > *Settings* > *Chamber configuration* (chap. 12.10).

7. Configuration of optional equipment

The "Optional equipment" menu item is visible only with optional chamber equipment.

To access the selection menu, go to Menu > Optional equipment

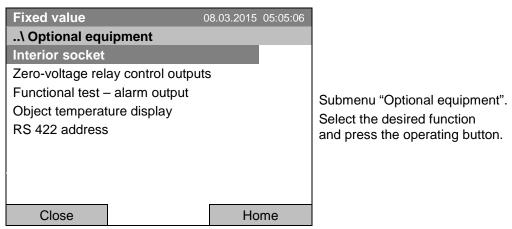

7.1 Turning on / off the optional interior socket

For chambers equipped with the water-protected interior socket (option, chap. 19.4) you can turn on and off the voltage of the interior socket via the controller.

To access the setting menu, go to Menu > Optional equipment > Interior socket

| Fixed value | 08.03.2015 05:05:06 | |
|----------------------|---------------------|---|
| \ Interior socket | | |
| Interior socket: Off | | Submenu "Interior socket". |
| | | The current setting is indicated. |
| | | To change the setting, press the operating button. |
| | | The modified setting is displayed. |
| | | "Interior socket: Off" = socket voltage-free "Interior socket: On" = socket with activated voltage |
| Close | Home | |
| | 1 | |

Go back to the initial view with "Home".

This symbol on the controller display indicates that the interior socket is activated.

7.2 Switching on or off the optional zero-voltage relay control outputs

For chambers equipped with zero-voltage relay outputs (available via BINDER INDIVIDUAL Customized Solutions, chap. 19.6), you can switch on or off the output via the controller.

To access the setting menu for the operating modes "Fixed value" and "Control off", go to Menu > Optional equipment > Zero-voltage relay control outputs.


| The setting for program op | eration is done through tr | le program eullor (chap. 9.3.7). | |
|----------------------------|----------------------------|---|--|
| Fixed value | 08.03.2015 05:05:06 | | |
| \ Zero-voltage relay co | ntrol outputs | Submenu "Zero-voltage relay control outputs". | |
| Control output 1: Off | | The current switching state of the zero-voltage rela | |
| Control output 2: Off | | control outputs is indicated. To change it, select the desired control output | |
| Control output 3: Off | | and press the operating button. | |
| | | The modified switching state is displayed. | |
| | | | |
| | | "Control output: Off" = zero-voltage relay output de- activated | |
| | | "Control output: On" = zero-voltage relay output acti- vated | |

The setting for program operation is done through the program editor (shap, 0.2.7)

Home

Go back to the initial view with "Home".

Close

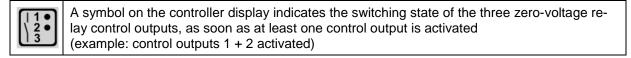

7.3 Functional test of the optional alarm output

For chambers equipped with the zero-voltage relay alarm output (option, chap. 19.3), you can switch on the output for test purpose via the controller and then switch it off again.

To access the setting menu, go to *Menu > Optional equipment > Functional test – alarm output*

| Fixed value | 08.03.2015 | 05:05:06 | | |
|----------------------------------|------------|----------|--|--|
| \ Functional test – alarm output | | | Submenu "Functional test – alarm output". | |
| Alarm output: Inactive | | | The current switching state is indicated. | |
| | | | To change the switching state, press the operating button. | |
| | | | The modified switching state is displayed. | |
| | | | "Alarm output: Inactive" = alarm output switched off "Alarm output: Active" = alarm output switched on (alarm state) | |
| Close | Hor | me | | |

Go back to the initial view with "Home".

7.4 Switching on or off the optional object temperature display

For chambers equipped with the digital object temperature display with a flexible Pt 100 temperature sensor (option, chap. 19.5), you can switch on or off the object temperature indication via the controller.

To access the setting menu, go to Menu > Optional equipment > Object temperature display

| Fixed value | 08.03.2015 05:05:06 | | |
|--------------------------------|---------------------|---|--|
| \ Object temperature display | | Submenu "Object temperature display". The current setting is indicated. To change the setting, press the operating button. | |
| object temperature display. On | | | |
| | | | |
| | | "Object temperature display: Off" = Object tempera- ture display deactivated | |
| | | "Object temperature display: On" = Object tempera- ture display activated | |
| Close | Home | | |

Go back to the initial view with "Home".

8. Set-point entry in "Fixed value" operating mode

8.1 Set-point ranges

| Temperature | Control and setting range: -10 °C / 14 °F up to +100 °C / 212 °F |
|-------------|--|
| Fan speed | 40 % up to 100 % (full speed). Reduce the fan speed only if required, because the spatial distribution of temperature will also be reduced. |
| | Technical data refers to 100% fan speed. |
| | |

| A | When you changed the temperature set-point, check the setting of the overtemperature safety controller class 3.1 (chap. 17.2) or the over-/undertemperature safety controller class 3.3 (option, chap. 17.3). |
|----------|---|
| | With set-point type "Limit", adapt the safety controller always when changing the temperature set-point. |

8.2 Entering the set-points via "quick menu"

| Fixed value \ Quick menu Measurement chart Active alarms Temperature setpoint Fan speed setpoint Safety controller setpoint Time second | 08.03.2015 05:05:06 | "Quick menu". Select the desired parameter and press the operating button. |
|--|---------------------|--|
| Time program Week program Close | Home | |
| When trying to enter | a setpoint in the | Controller off" operating mode, a notific: |

To enter set-points via quick menu, go to Quick menu.

When trying to enter a setpoint in the "Controller off" operating mode, a notification window shows "Controller mode is OFF!". Press the operating button to confirm with "Ok" and change the operating mode to "Fixed value" (chap. 6.2.1).

G

Temperature setting

To enter the temperature setpoint, go to Quick menu > Temperature setpoint

| Fixed value | 30 | 3.03.2015 05:05:06 | |
|---------------|----------|--------------------|--|
| \ Temperature | setpoint | | |
| 25.0 | 1 2 345 | [°C] | Entry n Select press t Setting Press t |
| Close | Ok | Home | |


Entry menu "Temperature setpoint". Select each number with the operating button and press the operating button to confirm. Setting range: -10 °C / 14 °F up to +100 °C / 212 °F Press the "Ok" button to confirm.

When entering a value outside the setting range, the message "invalid value" appears. Press the operating button to confirm with "Ok" and repeat the entry with a correct value.

Go back to the initial view with "Home" or enter the fan speed.

Fan speed setting

To enter the fan speed setpoint, go to Quick menu > Fan speed setpoint

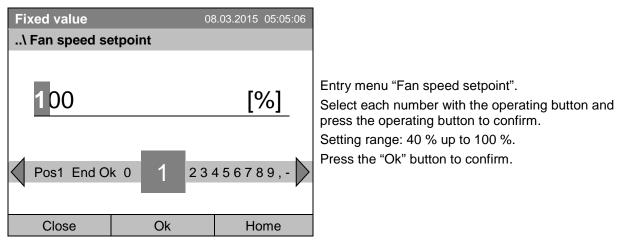

When entering a value outside the setting range, the message "invalid value" appears. Press the operating button to confirm with "Ok" and repeat the entry with a correct value.

Go back to the initial view with "Home".



When operating the fan with less than 100 % speed, the temperature performance and the spatial exactitude of the temperature may differ from the manufacturer's specifications. Do reduce the fan speed only if absolutely necessary due to special requirements.

8.3 Entering the set-points via general menu

To enter set-points via general menu, go to Menu > Setpoints

| Fixed value | 08.03.2015 | 05:05:06 | |
|-------------|------------|----------|---|
| \ Setpoints | | _ | |
| Temperature | | | |
| Fan speed | | | |
| | | | |
| | | | Submenu "Setpoints". |
| | | | Select "Temperature" or "Fan speed" and press the operating button. |
| | | | |
| | | | |
| | | | |
| Close | Ho | me | |

Temperature setting

To enter the temperature setpoint, go to *Menu* > *Setpoints* > *Temperature*

| Fixed value | | 08.03.2015 | 05:05:06 |
|---------------|----------|------------|-----------|
| \ Temperature | setpoint | | |
| 25.0 | | [°(| <u>C]</u> |
| Pos1 End Ok | 012 | 3456789 | 9,- |
| Close | Ok | Ho | me |

Entry menu "Temperature setpoint". Select each number with the operating button and press the operating button to confirm. Setting range: -10 °C / *14* °*F* up to +100 °C / *212* °*F* Press the "Ok" button to confirm.

When entering a value outside the setting range, the message "Invalid value" appears. Press the operating button to confirm with "Ok" and repeat the entry with a correct value.

Go back to the initial view with "Home" or enter the fan speed.


Fan speed setting

To enter the fan speed setpoint, go to Menu > Setpoints > Fan speed

| Fixed value | 30 | 3.03.2015 05:05:06 | |
|----------------|-----------|--------------------|---|
| \ Fan speed se | etpoint | | |
| 100 | k 0 1 234 | [%] 456789 | Entry menu "Fan speed setpoint". Select each number with the operating button and press the operating button to confirm. Setting range: 40 % up to 100 %. Press the "Ok" button to confirm. |
| Close | Ok | Home | |

When entering a value outside the setting range, the message "invalid value" appears. Press the operating button to confirm with "Ok" and repeat the entry with a correct value.

Go back to the initial view with "Home".

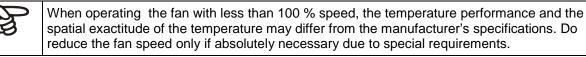

9. Time programs

The T4.12 program controller permits programming temperature cycles. It offers 52 program memory positions with up to 100 program sections each.

To access the menu selection for time programs, select Menu > Programs > Time program

| Fixed value | 0 | 8.03.2015 05:05:06 | |
|---------------------------|------------|--------------------|---|
| | Setpoint | Actual value | |
| Temperature [°(Fan [% | | | Initial view. Press the "Menu" button. |
| | | | |
| User | Quick menu | Menu | |




| Fixed value | 0 | 08.03.2015 05:05:06 | |
|--|-------------|---------------------|--|
| \ Menu | | | |
| Controller mode | | | |
| Event list | | | |
| Alarms | | | |
| | | | General menu. |
| Setpoints | | | Select "Programs" |
| Safety controller | r | | and press the operating button |
| Programs | | | and press the operating battern |
| Import/Export | | | |
| Settings | | $\overline{}$ | |
| 0 | | • | |
| Close | | Home | |
| 01030 | | Tiome | |
| Fixed value | 0 | 8.03.2015 05:05:06 | |
| \ Programs | | | |
| | _ | | |
| Time program | | | |
| Week program | | | |
| | | | Submenu "Programs". |
| | | | Select "Time Program" |
| | | | and press the operating button |
| | | | |
| | | | |
| | | | |
| | 1 | | |
| Close | | Home | |
| | | | |
| | | | |
| Fixed value | 0 | 08.03.2015 05:05:06 | |
| Fixed value \ Programs\Ti | | 08.03.2015 05:05:06 | |
| \ Programs\Ti | | 8.03.2015 05:05:06 | |
| \ Programs\Ti Start | | 08.03.2015 05:05:06 | |
| \ Programs\Ti Start Stop | | 08.03.2015 05:05:06 | |
| \ Programs\Tit Start Stop Pause | | 08.03.2015 05:05:06 | Submenu "Time programs". |
| \ Programs\Tin Start Stop Pause Resume | | 08.03.2015 05:05:06 | |
| \ Programs\Tit Start Stop Pause Resume Edit | | 08.03.2015 05:05:06 | |
| \ Programs\Tin Start Stop Pause Resume | | 08.03.2015 05:05:06 | Turn the operating button to see additional menu |
| \ Programs\Tit Start Stop Pause Resume Edit | | 08.03.2015 05:05:06 | Turn the operating button to see additional menu |
| \ Programs\Tit Start Stop Pause Resume Edit Create | | 08.03.2015 05:05:06 | Turn the operating button to see additional menu |
| \ Programs\Tit Start Stop Pause Resume Edit Create Rename | | 08.03.2015 05:05:06 | Turn the operating button to see additional menu |
| \ Programs\Tit Start Stop Pause Resume Edit Create Rename | | | Turn the operating button to see additional menu |
| \ Programs\Tit Start Stop Pause Resume Edit Create Rename Delete | | N8.03.2015 05:05:06 | Turn the operating button to see additional menu |
| \ Programs\Tit Start Stop Pause Resume Edit Create Rename Delete | me programs | | Turn the operating button to see additional menu |
| \ Programs\Tit Start Stop Pause Resume Edit Create Rename Delete Close Fixed value | me programs | Home | Turn the operating button to see additional menu |
| \ Programs\Tin Start Stop Pause Resume Edit Create Rename Delete Close Fixed value \ Programs\Tin | me programs | Home | Turn the operating button to see additional menu |
| \ Programs\Tit Start Stop Pause Resume Edit Create Rename Delete Close | me programs | Home | Turn the operating button to see additional menu |
| \ Programs\Tin Start Stop Pause Resume Edit Create Rename Delete Close Fixed value \ Programs\Tin | me programs | Home | Turn the operating button to see additional menu |
| \ Programs\Tin Start Stop Pause Resume Edit Create Rename Delete Close Fixed value \ Programs\Tin | me programs | Home | Turn the operating button to see additional menu |
| \ Programs\Tin Start Stop Pause Resume Edit Create Rename Delete Close Fixed value \ Programs\Tin | me programs | Home | Turn the operating button to see additional menu |
| \ Programs\Tin Start Stop Pause Resume Edit Create Rename Delete Close Fixed value \ Programs\Tin | me programs | Home | Turn the operating button to see additional menu items. |
| \ Programs\Tin Start Stop Pause Resume Edit Create Rename Delete Close Fixed value \ Programs\Tin | me programs | Home | Turn the operating button to see additional menu items. |
| \ Programs\Tin Start Stop Pause Resume Edit Create Rename Delete Close Fixed value \ Programs\Tin | me programs | Home | Turn the operating button to see additional menu items. |
| \ Programs\Tin Start Stop Pause Resume Edit Create Rename Delete Close Fixed value \ Programs\Tin | me programs | Home | Turn the operating button to see additional menu items. |
| \ Programs\Tin Start Stop Pause Resume Edit Create Rename Delete Close Fixed value \ Programs\Tin | me programs | Home | Turn the operating button to see additional menu items. |

9.1 Starting and running an existing time program

To start a time program, go to *Menu > Programs > Time program > Start*. (You can also go to *Quick menu > Time program > Start*, see below).

Starting is also possible directly from the program editor (chap. 9.3.9).

| (A) | In the "Control off" mode or during a running time or week program, no time program can be started. |
|-----|---|
| | |

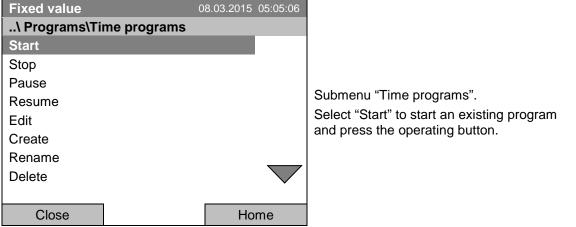

If no program has been created and saved so far, the message "No programs found" appears. Press the operating button to confirm with "Ok" and enter a program with "Create".

| Fixed value | 0 | 8.03.2015 05:0 | 5:06 | |
|-----------------|------------|----------------|------|--|
| \ Select progra | am | | | |
| Program0001 | | | | |
| Program0002 | | | | |
| Program0003 | | | | Submenu "Select program" (example). Select one of the programs and press the operating button to run the program |
| Close | | Home | | |
| Fixed value | 08 | 3.03.2015 05:0 | 5:06 | |
| \ Start date (D | D.MM.YYYY) | | | |
| 08.03.2 | | 3456789 | 9 | Entry menu "Start date". The current date is shown. For a postponed start, enter the desired start date with the operating but- ton. Press the "Ok" button to confirm. |
| Close | Ok | Home | | |



| Fixed value 08.03.2015 05:05:06 \ Start time (HH:MM:SS) | |
|--|--|
| 05:05:36 | Entry menu "Start time". The current time plus 30 seconds is shown. For a postponed start, enter the desired start time with the operating button. Press the "Ok" button to confirm. |
| Ins Pos1 End 0 1 2 3 4 5 6 7 8 9 Close Ok Home Time program 08.03.2015 05:05:06 | |
| Program0001 Setpoint Actual value Temperature [°C] 10.0 11.5 Section 20 End: 08.03.2015 08:07:44 100.0 Fan [%] 100.0 | Initial view (example values). The time program is running. |
| User Quick menu Menu | |



This symbol on the controller display indicates that a time program is running.

During a running time program, it is impossible to edit, rename or delete this program (when selecting such a function, a corresponding message is displayed). The other program functions are available.

During a running time program, no week program can be started.



To start a time program, you can also go to Quick menu > Time program > Start

| Fixed value | 08.03.2015 05:05:06 | |
|----------------------------|---------------------|---|
| \ Quick menu | | |
| Measurement chart | | |
| Active Alarms | | |
| Temperature setpoint | | |
| Fan speed setpoint | | "Quick menu". |
| Safety controller setpoint | | Select "Time program" |
| Time program | | and press the operating button |
| Week program | | |
| | | |
| | | |
| Close | Home | |
| | | 1 |
| Fixed value | 08.03.2015 05:05:06 | |
| \ Time program | | |
| | | |
| Start | | |
| | _ | |
| Start | _ | |
| Start | | Submenu "Time program". |
| Start | | Select "Start" to start an existing program |
| Start | | Submenu "Time program". Select "Start" to start an existing program and press the operating button. |
| Start | | Select "Start" to start an existing program |
| Start | | Select "Start" to start an existing program |
| Start | | Select "Start" to start an existing program |

The further procedure is equal to that described for the general menu.

Home

| Fixed value | 08.03.2015 | 05:05:06 | |
|------------------|------------|----------|---|
| \ Select program | | | |
| Program0001 | | | |
| Program0002 | | - | |
| Program0003 | | | |
| | | | Submenu "Select program" (example). |
| | | | Select one of the programs |
| | | | and press the operating button to start the program |
| | | | |
| | | | |
| | | | |
| Close | Но | me | |

Close

Performance after completing the program

The controller automatically changes to the "Fixed value" operation mode.

Before starting the program, check the temperature setpoint entered in the "Fixed value" operation mode. After end of the program, the temperature will equilibrate to this value.



NOTICE

Danger of damages by too high or too low temperature after the program ends. Damage to the charging material.

Before starting the program, check the temperature setpoint of "Fixed value" operation and if necessary adapt it.

9.2 Cancelling a running time program

To cancel a running time program, go to *Menu > Programs > Time program > Stop*.

To cancel a running time program, you can also go to Quick menu > Time program > Stop.

The controller returns to the initial view.

Performance after manual program stop

The controller automatically changes to the "Fixed value" operation mode.

Before starting a program, check the temperature setpoint entered in the "Fixed value" operation mode. After cancelling the program, the temperature will equilibrate to this value.



NOTICE

Danger of damages by too low temperature after cancelling the program. Damage to the charging material.

Before starting the program, check the temperature setpoint of "Fixed value" operation and if necessary adapt it.

9.3 Creating a new time program

For each program section you can enter a temperature set-point, the fan speed, the section's duration, the type of temperature transition "R" (ramp) or "S" (step) (see chap. 9.3.6), and the tolerance range.

| When changing the temperature set-point, check the setting of the safety controller (chap. 17.2 and 17.3) if the safety controller has been set to "Limit" mode. |
|--|
| |

If the fan is operated with less than 100 % speed, the temperature performance and the spatial exactitude of the temperature can differ from the manufacturer's specifications. The fan speed rate should only be reduced if absolutely necessary to meet special requirements.

Programming is saved in case of a power failure or after turning off the chamber.

To enter a new time program, go to *Menu > Programs > Time programs > Create*

| Fixed value | 08.03.2015 | 05:05:06 | |
|--------------------------|------------|---------------|---------------------------------|
| \ Programs\Time programs | | | |
| Start | | | |
| Stop | | | |
| Pause | | | • • • • • • |
| Resume | | | Submenu "Time program". |
| Edit | | | Select "Create" |
| Create | | | and press the operating button. |
| Rename | | | |
| Delete | | $\overline{}$ | |
| | | • | |
| Close | Ho | me | |

Creating a new program

| Fixed value | 08 | .03.2015 0 | 05:05:06 | |
|----------------|---------|------------|----------|--|
| \ Create new p | orogram | | | |
| New | | | | |
| Based on | | | | |
| | | | | Submenu "Create new program". |
| | | | | Select "New", to create an entirely new program, or "Based on…", to use an existing program for further editing and press the operating button. |
| Close | | Hom | ne | |

If you selected "Based on", then the program selection window appears:

| Fixed value \ Select program Program0001 Program0002 Program0003 | 08.03.2015 | 05:05:06 | |
|--|------------|----------|--|
| | | | Submenu "Select program" (example). Select the desired program and press the operating button. |
| Close | Ho | me | |

If no program has been created and saved so far, the message "No programs found" appears. Press the operating button to confirm with "Ok" and enter a program with "Create".

Now you can specify the name of the new time program:

| Enter name | 08 | 8.03.2015 05:05:06 | |
|------------|----------|--------------------|--|
| \ Program | | | |
| | | | |
| TUVWXY | (Z () 12 | 3456789 | Entry menu "Program". Enter the desired start time with the operating but- ton. Press the "Ok" button to confirm. |
| Close | Ok | Home | |

The time program editor is displayed. Following the selection "Based on...", this table for program entry shows the values of the selected program. Following the selection "New" there is an empty table, which can be filled section by section with sample values. You can edit the displayed values.

Entering the program values for the first parameter (temperature)

A first program line is shown. This corresponds to a program section. You can now edit the values.

| Fiz | xed va | lue | | | | 08.03.2015 05:05:06 | | | |
|-----|--------------------------|----------|------|------|---------|---------------------|-----------|-------|--|
| \ | \ Temperature controller | | | | | | | | |
| No. | Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S | 123 | |
| 1 | 25.00 | 00:00:15 | 1 | 0 | -999.00 | 999.00 | Ramp | 0 0 0 | |
| | | | | | | | | | Time program editor |
| | | | | | | | | | (view with optional control outputs). |
| | | | | | | | | | To edit the values, press the operating button, select |
| | | | | | | | | | the desired value and press the operating button |
| | | | | | | | | | again. |
| | | | | | | | | | |
| | | | | | | | | | |
| | ^ | | | L | | L | | | |
| | | | | Men | u | | \bigvee | | |

BINDER

To create a second program line (section), turn the operating button to the right and press it. The next section will be added.

| Fi | Fixed value 08 | | | | 08.03.2 | 015 05:0 | 05:06 | | |
|-----|--------------------------|----------|------|--------------------|---------|----------|-------|-----|--|
| \ | \ Temperature controller | | | | | | | | |
| No. | Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S | 123 | |
| 1 | 25.00 | 00:00:15 | 1 | 0 | -999.00 | 999.00 | Ramp | 000 | |
| 2 | 25.00 | 00:00:15 | 1 | 0 | -999.00 | 999.00 | Ramp | 000 | Time program editor |
| 3 | 25.00 | 00:00:15 | 1 | 0 | -999.00 | 999.00 | Ramp | 000 | (view with optional control outputs). |
| | | | | | | | | | To edit the values, press the operating button, select |
| | | | | | | | | | the desired value and press the operating button |
| | | | | | | | | | again. |
| | | | | | | | | | |
| | | | | | | | | | |
| | Menu | | | \bigtriangledown | | | | | |

The right column for the 3 control outputs is visible only on chambers equipped with optional control outputs.

9.3.1 Section handling

| Fixed value | | | Time program editor. | | | |
|------------------|-------|------|----------------------|----------|---------|---|
| \ Temperature co | ontro | ller | | | | |
| No. Value H:M:S | Ref. | Rep. | T. min | T.max | R/S | Select a value under "No." and press the operating button. |
| 1 25.00 00:00:15 | 1 | 0 | -999.00 | 999.00 | Ramp | and press the operating button. |
| \bigcup | | | | | | |
| Fixed value | | | 08.03 | .2015 05 | 5:05:06 | |
| \ Program sectio | ns | | | | | |
| Insert | | | | | | |
| Сору | | | | | | |
| Paste | | | | | | Submenu "Program sections". |
| Delete | | | | | | Select the desired function |
| | | | | | | and press the operating button. |
| | | | | | | |
| | | | | | | |
| | | | | | | |
| Close | | | | | | |

Note: When selecting "Insert" or "Paste", the new program section is inserted **before** the current section.

9.3.2 Temperature setpoint

| Fiz | ked value | | | 08.03.20′ | 15 05: | 05:06 | Time program editor. | | | | | |
|-----------------|--------------|---------|--------|-----------|--------|----------|---|--|--|--|--|--|
| \ | Temperature | contro | ller | | | | | | | | | |
| No. | Value H:M:S | S Ref. | Rep. T | min T. | max | R/S | Select a value under "Value" | | | | | |
| 1 | 25.00 00:00: | 15 1 | 0 -9 | 99.00 99 | 9.00 | Ramp | and press the operating button. | | | | | |
| | \bigcup | | | | | | | | | | | |
| Fiz | ked value | | | 08.03.20′ | 15 05: | 05:06 | | | | | | |
| \ | Temperature | setpoir | nt | | | | | | | | | |
| | 25.00 | | | [| °C] | <u> </u> | Entry menu "Temperature setpoint". A temperature value is shown. Enter the desired value with the operating button. | | | | | |
| \triangleleft | Pos1 End Ok | 01 | 2 3 | 4567 | 89, | - | Setting range: -10 °C / 14 °F up to +100 °C / 212 °F. Press the "Ok" button to confirm. | | | | | |
| | Close | (| Эk | | | | | | | | | |

When entering a value outside the setting range, the message "invalid value" appears. Press the operating button to confirm with "Ok" and repeat the entry with a correct value.

9.3.3 Section duration

| Fiz | Fixed value 08.03.2015 05:05:06 | | | | | | | Time program editor | | | |
|-----|---|--------------------|-------|------|---------|----------|---------|---|--|--|--|
| \ | Tempe | erature co | ntrol | ler | | | [°C] | Time program editor. Select a value under "H:M:S" | | | |
| No. | Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S | and press the operating button. | | | |
| 1 | 25.00 | 00:00:15 | 1 | 0 | -999.00 | 999.00 | Ramp | | | | |
| | \bigcup | | | | | | | | | | |
| Fiz | xed val | lue | | | 08.03 | .2015 05 | 5:05:06 | | | | |
| \ | Sectio | n duratio | n | | | | | | | | |
| 4 | Ins Pos | 00:15 s1 End Of | (| | 1234 | 5678 | 9 | Entry menu "Section duration". A duration value is shown. Enter the desired value with the operating button. Press the "Ok" button to confirm. | | | |
| | Close | e | (|)k | | | | | | | |

9.3.4 Repeating one or several sections within a time program

Enter the number of the target section, which shall be the start of the repeat cycle, under "Ref." and the number of repeats under "Rep.". To have sections repeated infinitely enter the number of cycles "Rep." as "-1".

| Fixed value 08.03.2015 05:05:06 \ Temperature controller [°C] No. Value H:M:S Ref. Rep. T. min T.max R/S 1 25.00 00:00:15 1 0 -999.00 999.00 Ramp | Time program editor. Select a value under "Ref." and press the operating button. | | |
|--|---|--|--|
| Fixed value 08.03.2015 05:05:06\ Reference section 1 | Entry menu "Reference section". Enter the target section number of the repeat cycle | | |
| Pos1 End Ok 1 2 3 4 5 6 7 8 9 , Close Ok | with the operating button. Press the "Ok" button to confirm. | | |
| Continue to enter the number of repeats: Fixed value 08.03.2015 05:05:06 Temperature controller [°C] No. Value H:M:S Ref. Rep. T. min T.max R/S 1 25.00 00:00:15 1 0 -999.00 999.00 Ramp | Time program editor. Select a value under "Rep." and press the operating button. | | |
| Fixed value 08.03.2015 05:05:06\ Number of repeats 0 | Entry menu "Number of repeats". Enter the desired number of repeats with the operat- | | |
| Del Pos1 End Ok 1 2 3 4 5 6 7 8 9 Close Ok | Press the "Ok" button to confirm. | | |


| Fiz | xed va | lue | | | 08.0 | 3.2015 0 | 5:05:06 | |
|-----|------------------|-----------|-------|-------|---------|----------|---------|----------------------|
| \ | Temp | erature o | contr | oller | | | [°C] | |
| No. | Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S | |
| 1 | 40.00 | 00:30:00 | 1 | 0 | -999.00 | 999.00 | Ramp | |
| 2 | 60.00 | 01:30:00 | 1 | 0 | -999.00 | 999.00 | Ramp | |
| 3 | 80.00 | 01:00:00 | 2 | 30 | -999.00 | 999.00 | Ramp | |
| 4 | 20.00 | 03:20:00 | 1 | 0 | 999.00 | 999.00 | Ramp | Time program editor. |
| | | | | | | | | |
| | | | | | | | | |
| - | | | | | | | | |
| | | | | | | | | |
| | | | | | | | | |
| | \bigtriangleup | | N | lenu | | \sim | 7 | |

The following example shows a time program where the sections 2 and 3 shall be repeated 30 times:

Sections 2 and 3 will be executed in total 31 times; only then will the program continue.

9.3.5 Tolerance range


You can specify a tolerance range for each program section with different values for the tolerance minimum and maximum. When the actual value exceeds the given threshold, the program is interrupted. This is indicated on the display, see chap. 9.4. When the actual temperature is situated again within the entered tolerance limits, the program automatically continues. Therefore, the duration of the program may be extended due to the programming of tolerances.

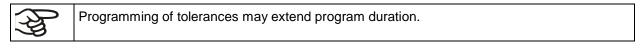

An entry of "-999" for the tolerance minimum means "minus infinite" and an entry of "999" for the tolerance maximum means "plus infinite". Entry of these values will never lead to program interruption.

When requesting a rapid temperature transition, we recommend not programming tolerance values in order to enable the maximum heating-up or cooling-down speed.

Start with the minimum value:

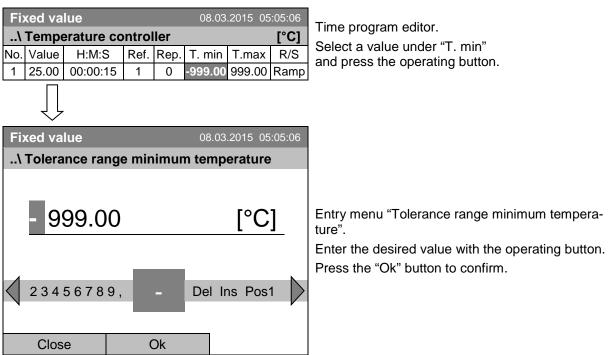

| F | Fixed value 08.03.2015 05:05:06 | | | | | | | Time program editor |
|----|---|---------------------|---------|------|---------|----------|--------|---|
| | Tempe | erature co | ontroll | er | | | [°C] | Time program editor. |
| No | . Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S | Select a value under "T. max" and press the operating button. |
| 1 | 25.00 | 00:00:15 | 1 | 0 - | 999.00 | 999.00 | Ramp | |
| | \int | | | | | | | _ |
| F | ixed va | lue | | | 08.03 | .2015 05 | :05:06 | |
| ' | Tolera | nce range | e max | imun | n temp | erature | e | |
| | 0123 | 9.00 3 4 5 6 7 8 | | | , - Del | [°C | | Entry menu "Tolerance range maximum tempera- ture". Enter the desired value with the operating button. Press the "Ok" button to confirm. |
| | Clos | е | 0 |)k | | | | |

Continue to enter the maximum value:

9.3.6 Set-point ramp and set-point step modes

"Ramp" mode

The set-point of a given program section functions as the section's target temperature. During the section's duration, the temperature set-point gradually passes from the set-point of the previous program section to this target set-point. The actual temperature value follows the continually changing set-point.

Programming in the "ramp" mode allows various kinds of temperature transitions:

Gradual slow temperature changes

The set-point changes its value gradually during the entered section duration. The actual temperature value (X) follows the continually moving set-point (W) at any time.

• Constant temperature

The initial values of two subsequent program sections are identical; so the temperature remains constant during the duration of the first program section.


Sudden temperature changes

These changes occur rapidly within the minimum amount of time (minimum entry: 1 second).

"Step" mode


The set-point of a given program section functions as the section's target temperature. At the start of the program section, the chamber heats up or cools down with the maximum speed to reach the entered value; and then it holds it for the remaining section time. The set-point temperature remains constant for the section's duration.

Selecting the setting "Ramp" or "Step":

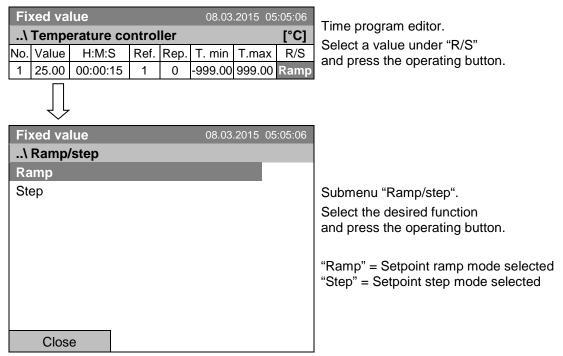

BINDER


Examples:

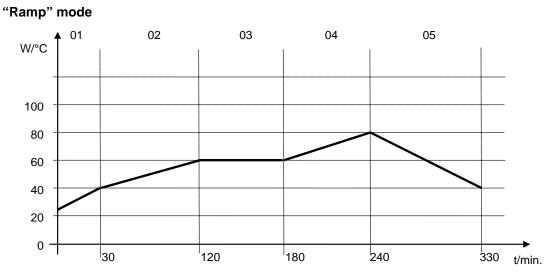

Corresponding program table:

| No. | Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S |
|-----|-------|----------|------|------|--------|-------|------|
| 01 | 40.0 | 00:30:00 | 1 | 0 | -999 | +999 | Ramp |
| 02 | 60.0 | 01:30:00 | 1 | 0 | -5 | +5 | Ramp |
| 03 | 60.0 | 01:00:00 | 1 | 0 | -2 | +2 | Ramp |
| 04 | 80.0 | 01:00:00 | 1 | 0 | -999 | +999 | Ramp |
| 05 | 40.0 | 01:30:00 | 1 | 0 | -999 | +999 | Ramp |

"Step" mode

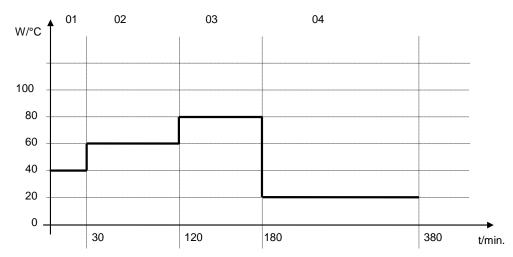

Corresponding program table

| No. | Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S |
|-----|-------|----------|------|------|--------|-------|------|
| 01 | 40.0 | 00:30:00 | 1 | 0 | -999 | +999 | Step |
| 02 | 60.0 | 01:30:00 | 1 | 0 | -5 | +5 | Step |
| 03 | 80.0 | 01:00:00 | 1 | 0 | -2 | +2 | Step |
| 04 | 20.0 | 03:20:00 | 1 | 0 | -999 | +999 | Step |

9.3.7 Switching on or off the optional zero-voltage relay outputs

For chambers equipped with zero-voltage relay outputs (option, chap. 19.6), you can switch on or off the outputs for each program section via the program editor.

| Fixed value 08.03.2015 05:05: \ Temperature controller [°(C) No. Value H:M:S Ref. Rep. T. min T.max R/S T 1 25.00 00:00:15 1 0 -999.00 999.00 Rampe I | Time program editor (with option relay outputs) Select a field under "123" and press the operating button |
|--|--|
| Fixed value 08.03.2015 05:05:06 Zero -voltage relay outputs Outputs 1 [Off] 2[Off] 3[Off] Outputs 1 [Off] 2[Off] 3[Off] Outputs 1[Off] 2[On] 3[Off] Outputs 1[Off] 2[On] 3[Off] Outputs 1[Off] 2[On] 3[Off] Outputs 1[Off] 2[On] 3[Off] Outputs 1[Off] 2[Off] 3[On] Outputs 1[Off] 2[Off] 3[On] Outputs 1[Off] 2[On] 3[On] Outputs 1[On] 2[On] 3[On] Outputs 1[On] 2[On] 3[On] Close | Submenu "Zero -voltage relay outputs". The possible combination of switching states are in- dicated. Select the desired combination of switching states and press the operating button." [On] = zero-voltage relay outputs turned on [Off] = zero-voltage relay outputs turned off |

 $\begin{bmatrix} 1 & \bullet \\ 1 & 2 & \bullet \\ 3 & 0 & 1 \\ 1$

A symbol on the controller display indicates the switching state of the three zero-voltage relay outputs as soon as at least one output is switched on (example: outputs 1 + 2 turned on)

9.3.8 Calling up the next parameter

| Fixed value 08.03.2 | | | | | | | 5:05:06 | |
|---------------------|-------|----------|-------|-------|---------|--------|---------|------------------------------|
| \ | Tempe | erature | contr | oller | | | [°C] | |
| No. | Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S | |
| 1 | 40.00 | 00:30:00 | 1 | 0 | -999.00 | 999.00 | Step | |
| 2 | 60.00 | 01:30:00 | 1 | 0 | -5.00 | 5.00 | Step | |
| 3 | 80.00 | 01:00:00 | 1 | 0 | -5.00 | 5.00 | Step | Time program editor (example |
| 4 | 20.00 | 03:20:00 | 1 | 0 | -999.00 | 999.00 | Step | Press the "Menu" button. |
| 5 | 40.00 | 00:30:00 | 1 | 0 | -999.00 | 999.00 | Step | |
| 6 | 60.00 | 01:30:00 | 1 | 0 | -5.00 | 5.00 | Step | |
| 7 | 70.00 | 01:00:00 | 1 | 0 | -5.00 | 5.00 | Step | |
| 8 | 60.00 | 01:00:00 | 1 | 0 | -5.00 | 5.00 | Step | |
| | | | Menu | | | | | |



| Fixed value | 08.03.2015 | 05:05:06 | |
|------------------------|------------|----------|---|
| \ Program menu | | | |
| Select parameter | | | |
| Save and exit | | | |
| Save and run | | | |
| Exit (without saving!) | | | Submenu "Program menu". Select "Select parameter" and press the operating button. |
| | | | |
| Close | | | |
| | | | |

(Ag

With the "Close" button, the controller returns to the time program editor.

| Fixed value | | 08.03.2015 | 05:05:06 | |
|---------------------|--------|------------|----------|------------------------------------|
| \ Program param | neters | | | |
| Temperature control | oller | | | |
| Fan speed | | | | |
| | | | | Submenu Select "Fa and press |
| Close | | | | |

Submenu "Program parameters". Select "Fan speed main controller" and press the operating button.

Entering the set-point values for another parameter (fan speed)

The number of program lines (program sections) equal to the number in the temperature program is displayed. The settings of section length, repeats and the selection "Ramp" or "Step" are taken over from the temperature program; they are no editable in this view. The symbol " *** " indicates that there are no tolerance value required for the fan speed. You can enter the fan speed set-points.

If you want to insert further sections or perform any other programming, first change back to the temperature program through *Menu* > *Select parameter* > *Temperature controller*.

| Fi | xed va | lue | | | 08.0 | 3.2015 0 | 5:05:06 |
|-----|------------------|----------|------|------|--------|---------------------|---------|
| \ | Fan sp | beed | | | | | [%] |
| No. | Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S |
| 1 | 100.00 | 00:30:00 | 1 | 0 | *** | *** | Ramp |
| 1 | 100.00 | 00:10:00 | 1 | 0 | *** | *** | Ramp |
| 1 | 100.00 | 00:30:00 | 2 | 3 | *** | *** | Ramp |
| 1 | 100.00 | 01:00:00 | 1 | 0 | *** | *** | Ramp |
| | | | | | | | |
| | | | | | | | |
| | | | | | | | |
| | | | | | | | |
| | | | | | | | |
| | \bigtriangleup | | Μ | lenu | | $\overline{\nabla}$ | 7 |

Entering the fan speed setpoint

| Fixed value | | 08.0 | 3.2015 0 | 5:05:06 | |
|-------------------|-----------|--------|----------|-----------|---|
| \ Fan speed [%] | | | | | Time program editor. |
| No. Value H:M:S | Ref. Rep. | T. min | T.max | R/S | Select a value under "Value" and press the operating button. |
| 1 100.00 00:00:15 | 1 0 | *** | *** | Ramp | |
| \bigcup | | | | | |
| Fixed value | | 08.0 | 3.2015 0 | 5:05:06 | |
| \ Fan setpoint | | | | | |
| 100 | _ | _ | [% | <u>5]</u> | Entry menu "Fan setpoint". A fan speed value is shown. Enter the desired value with the operating button. Setting range: 40 % up to 100 % Press the "Ok" button to confirm. |
| Pos1 End Ok | 0 1 Ok | 234 | 5678 | 9, | |

When entering a value outside the setting range, the message "invalid value" appears. Press the operating button to confirm with "Ok" and repeat the entry with a correct value.

9.3.9 Saving the time program and leaving the program editor

| Fiz | xed va | lue | | | 5:05:06 | | | |
|-----|-------------------------------|-----------|------|------|---------|----------|---------------------------------|--------------------------------|
| \ | \ Temperature controller [°C] | | | | | | | |
| No. | Value | H:M:S | Ref. | Rep. | T. min | T.max | R/S | |
| 1 | 40.00 | 00:30:00 | 1 | 0 | -999.00 | 999.00 | Step | |
| 2 | 60.00 | 01:30:00 | 1 | 0 | -5.00 | 5.00 | Step | |
| 3 | 80.00 | 01:00:00 | 1 | 0 | -5.00 | 5.00 | Step | Time program editor (example). |
| 4 | 20.00 | 03:20:00 | 1 | 0 | -999.00 | 999.00 | Step | Press the "Menu" button. |
| 5 | 40.00 | 00:30:00 | 1 | 0 | -999.00 | 999.00 | Step | |
| 6 | 60.00 | 01:30:00 | 1 | 0 | -5.00 | 5.00 | Step | |
| 7 | 70.00 | 01:00:00 | 1 | 0 | -5.00 | 5.00 | Step | |
| 8 | 60.00 | 01:00:00 | 1 | 0 | -5.00 | 5.00 | Step | |
| | \bigtriangleup | | N | lenu | | \sim | 7 | |
| | | | | | | | | 1 |
| Fiz | xed va | lue | | | 08.0 | 3.2015 0 | 5:05:06 | |
| \ | Progra | am meni | u | | | | | |
| Se | elect pa | arameter | | | | | | |
| Sa | ive and | d exit | | | | | | |
| Sa | ve and | d run | | | | | | |
| Fx | it (with | out savir | na!) | | | | | Submenu "Program menu". |
| | | | ·9·/ | | | | | Select the desired function |
| | | | | | | | and press the operating button. | |
| | | | | | | | | |
| | | | | | | | | |
| | | | | | | | | |
| | | | | | | | | |
| | Clos | е | | | | | | |



With the "Close" button, the controller returns to the time program editor.

| Select parameter | Changing between temperature and fan speed. | | |
|------------------------|--|--|--|
| Save and exit | Saving the program. The controller returns to the "Time programs" submenu. You can now select and start the program as described in chap. 9.1. | | |
| Save and run | Saving the program and start it. Enter the start date and time, see chap. 9.1. If another time or week program is running, the program is only saved, but not started. A corresponding message is displayed. | | |
| Exit (without saving!) | Attention: the program is not saved. After a security question, the controller returns to the initial view. | | |
| Button "Close" | The controller returns to the time program editor. You can continue programming. | | |

Make sure that you saved the time program before leaving the program editor.

With "Exit (without saving!)" you exit the program editor without saving the program. There is a security question first:

| Fixed value | 08.03.2015 | 05:05:06 | |
|-------------------------------|------------|----------|-------------|
| \ Exit (without saving) | | | |
| Do not exit | | | |
| Really exit (without saving!) | | | |
| | | | |
| | | | Submenu " |
| | | | This a secu |
| | | | and press t |
| | | | |
| | | | |
| | | | |
| Close | | | |
| | | | i |

Submenu "Exit (without saving)". This a security question. Select the desired function and press the operating button.

If you selected "Really exit (without saving!)", the controller goes back to the initial view.

9.4 **Program interruption**

You can manually interrupt a time program (pause), or this will automatically occur when exceeding the entered tolerance range values of the corresponding program section (see chap. 9.3.5).

Manual program interruption

To interrupt a time program, go to *Menu > Programs > Time programs > Pause*



| Time program | 08 | .03.2015 05:05:06 | |
|-------------------|----------------------|----------------------|--|
| \ Time program | n | | |
| Start | | | |
| Stop | | | |
| Pause | | | Submenu "Time program". |
| Resume | | | Select "Pause" to interrupt the running time program |
| Edit | | | and press the operating button. |
| Create | | | |
| Rename | | | |
| Delete | | | |
| Close | | Home | |
| With button "Home | e" the controller re | eturns to the initia | al view. |
| Time program p | baused 08 | .03.2015 05:05:06 | |
| Program0001 | Setpoint | Actual value | |
| Temperature [°C | 10.0 | 11.5 | |

This symbol on the controller display indicates that a running time program is interrupted.

Continuing the time program after a manual interruption

| Time program | paused | 08.03.2015 | 05:05:06 |
|----------------|--------|------------|--------------------|
| \ Time program | n | | |
| Start | | | |
| Stop | | | |
| Pause | | | |
| Resume | | | |
| Edit | | | |
| Create | | | |
| Rename | | | |
| Delete | | | \bigtriangledown |
| | | | |
| Close | | Но | me |

Submenu "Time program". Select "Resume" to continue the time program and press the operating button.

With button "Home" the controller returns to the initial view.

9.5 Deleting a time program

To delete a time program, go to *Menu > Programs > Time program > Delete* or *Menu > Programs > Time program > Delete all*

| Fixed value | 0 | 8.03.2015 05:05:06 | |
|----------------|------------|--------------------|---|
| \ Programs\Tim | ne program | | |
| Start | | | |
| Stop | | | |
| Pause | | | |
| Resume | | | Submenu "Time program". |
| Edit | | | Select "Delete" or "Delete all" (next page) and press the operating button |
| Create | | | and press the operating button |
| Rename | | | |
| Delete | | | |
| | | | |
| Close | | Home | |

If you selected "Delete", select then the time program to be deleted and press the operating button.

If you selected "Delete all", all time programs will be deleted in the controller.

Before deleting there is a security question:

| Fixed value | C | 8.03.2015 | 05:05:06 | |
|----------------|--------|-----------|----------|--|
| \ Confirm oper | ration | | | |
| Do not delete | | | | |
| Really delete | | | | |
| | | | | Submenu "Confirm operation". This a security question. Select the desired function and press the operating button. |
| Close | | Hor | ne | |

Go back to the initial view with "Home".



10. Week programs

The T4.12 program controller permits programming week programs with real-time reference. It offers 8 week program places in total with up to 30 shift points for each week program.

To access the menu selection for week programs, select Menu > Programs > Week program

| Fixed value | 08 | 8.03.2015 05:05:06 | |
|-------------------|-------------|--------------------|---|
| | Setpoint | Actual value | |
| Temperature [°C | 25.0 | 25.2 | |
| Fan [%] | 100.0 | | Initial view. |
| | | | Press the "Menu" button. |
| | | | |
| ▶●◀ | | | |
| User | Quick menu | Menu | |
| Fixed value | 0 | 3.03.2015 05:05:06 | |
| \ Menu | | | |
| Controller mode | | | |
| Event list | | | |
| Alarms | | | General menu. |
| Setpoints | | | |
| Safety controller | | | Select "Programs" and press the operating button |
| Programs | | | |
| Import/Export | | | |
| Settings | | | |
| Close | | Home | |
| Fixed value | 08 | 3.03.2015 05:05:06 | |
| \ Programs | | | |
| Time program | | | |
| Week program | | | |
| | | | Submenu "Programs". |
| | | | Select "Week Program" |
| | | | and press the operating button |
| | | | |
| | | | |
| | | | |
| Close | | Home | |

10.1 Starting and running an existing week program

To start a week program, go to *Menu > Programs > Week program > Start*

(You can also go to Quick menu > Week program > Start, see below).

Starting is also possible directly from the program editor (chap. 10.3.8).

| A | In the "Control off" mostarted. | ode or during a runr | ning time or week program, no week program can be |
|--|--------------------------------------|----------------------|--|
| Fixed va \ Progra Start Stop Edit Create Rename Delete Delete al | ams\Week Programs | 08.03.2015 05:05:06 | Submenu "Week Programs". Select "Start" to start an existing week program and press the operating button |
| Fixed va \ Select Program Day-Nigh | program 0004 0005 It | 08.03.2015 05:05:06 | Submenu "Select program" (example). Select one of the programs and press the operating button to run the program |

If no program has been created and saved so far, the message "No programs found" appears. Press the operating button to confirm with "Ok" and enter a program with "Create".

| Fixed value | 30 | 3.03.2015 05:05:06 | |
|-----------------|------------|--------------------|--|
| \ Start date (D | D.MM.YYYY) | | |
| 0 8.03.2 | _ | 7 8 9 , - Del | Entry menu "Start date". The current date is shown. For a postponed start, enter the desired start date with the operating but- ton. Press the "Ok" button to confirm. |
| Close | Ok | Home | |



| Fixed value 08.03.2015 05:05:06 \ Start time (HH:MM:SS) | |
|---|--|
| 05:05:36 | Entry menu "Start time". The current time plus 30 seconds is shown. For a postponed start, enter the desired start time with the operating button. Press the "Ok" button to confirm. |
| Close Ok Home | |
| Week program08.03.201505:05:06Program0004SetpointActual value | |
| Temperature [°C] 15.0 15.3 Section: 0002 End: Sunday 00:00:00 | Initial view. |
| Fan [%] 100.0 | The week program is running. |
| User Quick menu Menu | |

This symbol on the controller display indicates that a week program is running.

During a running week program, it is impossible to edit, rename or delete this program (when selecting such a function, a corresponding message is displayed). The other program functions are available.

During a running week program, no time program can be started.

To start a week program, you can also go to Quick menu > Week program > Start

| Fixed value | 08.03.2015 | 05:05:06 | |
|----------------------------|------------|----------|--------------------------------|
| \ Quick menu | | | |
| Measurement chart | | | |
| Active Alarms | | | |
| Temperature setpoint | | | |
| Fan speed setpoint | | | "Quick menu". |
| Safety controller setpoint | | | Select "Week program" |
| Time program | | | and press the operating button |
| Week program | | | |
| | | | |
| | | | |
| | | | |
| Close | Но | me | |



| Fixed value | 08.03 | 3.2015 05:05:06 | |
|----------------|-------|-----------------|---|
| \ Week progran | n | | |
| Start | | | |
| Stop | | | |
| | | | |
| | | | Submenu "Week program". |
| | | | Select "Start" to start an existing program |
| | | | and press the operating button. |
| | | | |
| | | | |
| | | | |
| Class | | Homo | |
| Close | | Home | |

The further procedure is equal to that described for the general menu.

| Fixed value \ Select program Program0001 Program0002 Program0003 | 08.03.2015 05:05:06 | Submenu "Select program" (example). Select one of the programs and press the operating button to start the program |
|--|---------------------|--|
| Close | Home | |

10.2 Cancelling a running week program

To cancel a week program, go to *Menu > Programs > Week program > Stop*.


To cancel a running week program, you can also go to Quick menu > Week program > Stop.

The controller returns to the initial view.

Performance after manual program stop

The controller automatically changes to the "Fixed value" operation mode.

Before starting a program, check the temperature setpoint entered in the "Fixed value" operation mode. After cancelling the program, the temperature will equilibrate to this value.

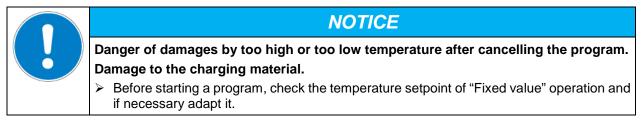

10.3 Creating a new week program

A week program permits defining up to 30 sections for the whole week. These sections function as shift points. A shift-point is characterized by its time, temperature value, and state (active / inactive). With a running week program, the temperature of the currently active shift point is maintained until the moment of the next active shift point with its new temperature set-point.

Example:

| No. | Value | Day | H:M:S | Activity |
|-----|-------|---------|----------|----------|
| 1 | 35.00 | Monday | 08:00:00 | Active |
| 2 | 40.00 | Monday | 10:00:00 | Active |
| 3 | 35.00 | Monday | 18:00:00 | Active |
| 4 | 10.00 | Monday | 20:00:00 | Active |
| 5 | 35.00 | Tuesday | 08:00:00 | Active |
| | | | | |

Programming is saved in case of a power failure or after turning off the chamber.

To enter a new week program, go to *Menu > Programs > Week program > Create*

| Fixed value\ Programs\Week programs Start Stop Edit Create Rename Delete Delete all | 08.03.2015 05:05:06 | Submenu "Week program". Select "Create" and press the operating button. |
|---|---------------------|--|
| Close | Home | |
| Fixed value Create new program New Based on | 08.03.2015 05:05:06 | Submenu "Create new program". Select "New", to create an entirely new program, or "Based on", to use an existing program for further editing and press the operating button. |
| Close | Home | |

If you selected "Based on", then the program selection window appears:

| Fixed value | 08.03.2015 | 05:05:06 | |
|------------------|------------|----------|---|
| \ Select program | | | |
| Program0004 | | | |
| Program0005 | | | |
| Day-Night | | | |
| | | | Submenu "Select program" (example). |
| | | | Select the desired program and press the operating button. |
| | | | and press the operating button. |
| | | | |
| | | | |
| | | | |
| Close | Ho | me | |

If no program has been created and saved so far, the message "No programs found" appears. Press the operating button to confirm with "Ok" and enter a program with "Create".

Now you can specify the name of the new week program:

| Enter name | 08 | 8.03.2015 05:05:06 | |
|------------|----------|--------------------|--|
| \ Program | | | |
| TUVWXY | ′Z () 12 | 3456789 | Entry menu "Program". Enter the desired start time with the operating but- ton. Press the "Ok" button to confirm. |
| Close | Ok | Home | |

The week program editor is displayed. Following the selection "Based on...", this table for program entry shows the values of the selected program. Following the selection "New" there is an empty table, which can be filled section by section with sample values. You can then edit the displayed values.

Entering the program values for the first parameter (temperature)

A first program line is shown. This corresponds to a program section. You can now edit the values.

| Fixe | d value | | 08. | 03.2015 05: | 05:06 | |
|------|-------------------------------|--------|----------|---------------------|-------|--|
| \ Te | \ Temperature controller [°C] | | | | | |
| No. | Value | Day | H:M:S | Activity | 123 | |
| 1 | 25.00 | Monday | 00:00:00 | Inactive | 0 0 0 | |
| | | | | | | Week program editor |
| | | | | | | (view with optional control outputs) |
| | | | | | | To edit the values, select the desired value and |
| | | | | | | press the operating button. |
| | | | | | | |
| | | | | | | |
| | | | | | | |
| | | Ме | nu | $\overline{\nabla}$ | | |

To create a second program line (section), turn the operating button to the right and press it. The next section will be added.

| Fixe | Fixed value 08.03.2015 05:05:06 | | | | | | | |
|------------------------------|---------------------------------|--------|----------|----------|-----|--|--|--|
| \ Temperature controller [°C | | | | | | | | |
| No. | Value | Day | H:M:S | Activity | 123 | | | |
| 1 | 25.00 | Monday | 00:00:00 | Inactive | 000 | | | |
| 2 | 25.00 | Monday | 00:00:00 | Inactive | 000 | | | |
| | | | | | | | | |
| | | | | | | | | |
| | | | | | | | | |
| | | | | | | | | |
| | | | | | | | | |
| | | | | | | | | |
| | \bigtriangleup | Ме | nu | \sim | | | | |


Week program editor (view with optional control outputs)

To edit the values, press the operating button, select the desired value and press the operating button again.

10.3.1 Section handling

| Fixe | d value | | 08.03.2 | 015 05:05:06 | | |
|--------|---------------------|--------------|----------------------|--------------|---|--|
| \ Te | mperatu | re controlle | Week program editor. | | | |
| No. | Value | Day | H:M:S | Activity | Select a value under "No." and press the operating button. | |
| 1 | 25.00 | Monday | 00:00:00 | Inactive | | |
| _ | $\bigcup_{i=1}^{n}$ | | | | | |
| Fixe | d value | | 08.03.2 | 015 05:05:06 | | |
| \ Pr | ogram se | ections | | | | |
| Inse | rt | | | | | |
| Сору | / | | | | | |
| Paste | e | | | | | |
| Delete | | | | | Submenu "Program sections". | |
| | | | | | Select the desired function | |
| | | | | | and press the operating button. | |
| | | | | | | |
| | | | | | | |
| | | | | | | |
| (| Close | | | | | |


Note: When selecting "Insert" or "Paste", the new program section is inserted **before** the current section.

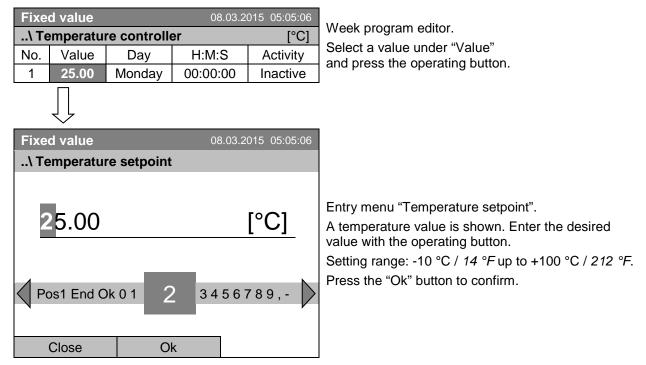

10.3.2 Temperature setpoint

When entering a value outside the setting range, the message "invalid value" appears. Press the operating button to confirm with "Ok" and repeat the entry with a correct value.

10.3.3 Day of the week

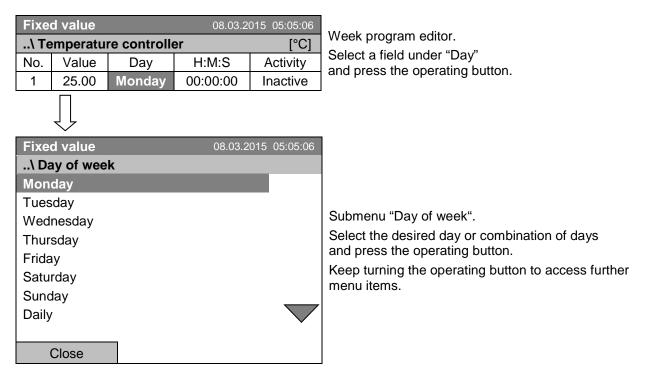

| Fixed value | 08.03.2015 05:05:06 | |
|---|---------------------|--|
| \ Day of week | | |
| Mon-Fri Sun-Thu Mon-Sat Fri+Sat Sat+Sun | | Submenu "Settings" (next page) Select the desired function and press the operating button. |
| Close | | |

10.3.4 Time of the day

| Fixed value 08.03.2015 05:05:0 | | | | 015 05:05:06 | | | |
|--------------------------------|----------|-------------|----------|--------------|---|--|--|
| \ Tem | nperatur | e controlle | er | [°C] | Week program editor. | | |
| No. | Value | Day | H:M:S | Activity | Select a value under "H:M:S" and press the operating button. | | |
| 1 | 25.00 | Monday | 00:00:00 | Inactive | and press the operating button. | | |
| | Ţ | | | | | | |
| Fixed | value | | 08.03.20 | 015 05:05:06 | | | |
| \ Tim | e of day | 1 | | | | | |
| Ins | Pos1 Er | | | 6789 | Entry menu "Time of day". Enter the desired moment of the shift-point with the operating button. Press the "Ok" button to confirm. | | |

| Fixe | d value | | | | |
|-------|---------------------|--------------|----------|--------------|---|
| \ Te | emperatu | re controlle | er | [°C] | Week program editor. |
| No. | Value | Day | H:M:S | Activity | Select a field under "Activity" and press the operating button. |
| 1 | 25.00 | Monday | 00:00:00 | Inactive | and press the operating button. |
| | $\bigcup_{i=1}^{n}$ | | | | |
| Fixe | d value | | 08.03.2 | 015 05:05:06 | |
| \ Ac | ctivity | | | | |
| Inact | tive | | | | |
| Activ | e | | | | |
| | | | | | Submenu "Activity". Select the desired function and press the operating button. |
| | Close | | | | |

10.3.5 Activity of the shift-point

10.3.6 Switching on or off the optional zero-voltage relay outputs

| Fixed value 08.03.2015 05:05:06 | | | | | | | | |
|---------------------------------|-------|--------|----------|----------|-----|----|--|--|
| \ Temperature controller [°C] | | | | | | | | |
| No. | Value | Day | H:M:S | Activity | 123 | S | | |
| 1 | 25.00 | Monday | 00:00:00 | Inactive | 000 | ar | | |
| | | | | | | | | |

Week program editor (view with optional control outputs) Select a field under "1123" and press the operating button.

| Fixed value | 08.03.2015 05:05:06 | | | | | | | | |
|----------------------------|-------------------------------|--|--|--|--|--|--|--|--|
| \ Zero -voltage | \ Zero -voltage relay outputs | | | | | | | | |
| Outputs 1 [Off] | 2[Off] 3[Off] | | | | | | | | |
| Outputs 1[On] 2 | [Off] 3[Off] | | | | | | | | |
| Outputs 1[Off] 2 | [On] 3[Off] | | | | | | | | |
| Outputs 1[On] 2 | [On] 3[Off] | | | | | | | | |
| Outputs 1[Off] 2 | [Off] 3[On] | | | | | | | | |
| Outputs 1[On] 2 | [Off] 3[On] | | | | | | | | |
| Outputs 1[Off] 2[On] 3[On] | | | | | | | | | |
| Outputs 1[On] 2[On] 3[On] | | | | | | | | | |
| | | | | | | | | | |
| Close | | | | | | | | | |

Submenu "Zero -voltage relay outputs".

The possible combination of switching states are indicated.

Select the desired combination of switching states and press the operating button."

[On] = zero-voltage relay outputs turned on [Off] = zero-voltage relay outputs turned off

 \downarrow

A symbol on the controller display indicates the switching state of the three zero-voltage relay outputs as soon as at least one output is switched on (example: outputs 1 + 2 turned on)

| Fixe | d value | | | | | |
|--------------------------|------------------|---------|------------------------|----|--------------------|--------------------------------|
| \ Temperature controller | | | | | | |
| No. | Value | Day | H:M:S | | Activity | |
| 1 | 35.00 | Monday | 08:00: | 00 | Active | |
| 2 | 40.00 | Monday | 10:00: | 00 | Active | |
| 3 | 35.00 | Monday | 18:00: | 00 | Active | Week program editor (example). |
| 4 | 10.00 | Monday | 20:00: | 00 | Active | Press the "Menu" button. |
| 5 | 35.00 | Tuesday | 08:00:00 | | Active | |
| 6 | 40.00 | Tuesday | 10:00: | 00 | Active | |
| 7 | 35.00 | Tuesday | 18:00: | 00 | Active | |
| 8 | 10.00 | Tuesday | 20:00: | 00 | Active | |
| | \bigtriangleup | Men | u | | \bigtriangledown | |
| | | | | | | 1 |
| Fixe | d value | | | | | |
| \ Pr | ogram m | enu | | | | |
| Sele | ct param | eter | | | | |
| Save | and exit | | | | | |
| Save | and run | | Submenu "Program menu" | | | |

10.3.7 Calling up the next parameter

Submenu "Program menu". Select "Select parameter" and press the operating button.

Close

Exit (without saving!)

With the "Close" button, The controller returns to the week program editor.

| | | | - |
|------------------------|------------|----------|-----------------------------|
| Fixed value | 08.03.2015 | 05:05:06 | |
| \ Program parameters | | | |
| Temperature controller | | | |
| Fan speed | | | |
| | | | Submonu "F |
| | | | Submenu "P |
| | | | Select "Fan and press th |
| | | | |
| | | | |
| | | | |
| | | | |
| Close | | | |

Entering the set-point values for another parameter (fan speed)

The number of program lines (program sections) equal to the number in the temperature program is displayed. The settings of the shift points (day, time, activity) are taken over from the temperature program; they are no editable in this view. You can enter the fan speed set-points.



If you want to insert further sections or perform any other programming, first change back to the temperature program through *Menu* > *Select parameter* > *Temperature controller*.

| Fixe | d value | | 08.03 | 3.2015 05:05:06 | |
|------|----------|--------|----------|--------------------|-------------------------------|
| \ Fa | an speed | | | [%] | |
| No. | Value | Day | H:M:S | Activity | |
| 1 | 100.00 | Monday | 10:00:00 |) Active | |
| 1 | 100.00 | Monday | 12:00:00 |) Active | |
| 1 | 100.00 | Monday | 00:00:00 |) Inactive | |
| | | | | | Week program editor (example) |
| | | | | | |
| | | | | | |
| | | | | |] |
| | | | | | |
| Menu | | | | \bigtriangledown |] |

Entering the fan speed setpoint

| Fixe | d value | | 08.03.201 | 5 05:05:06 | Week program editor |
|------|---------------|--------|-----------|---------------------|---|
| \ Fa | an speed | | | [%] | Week program editor. |
| No. | Value | Day | H:M:S | Activity | Select a value under "Value" and press the operating button. |
| 1 | 100.00 | Monday | 00:00:00 | Inactive | |
| | \bigcup | | | | |
| Fixe | d value | | 08.03.201 | 15 05:05:06 | |
| \ Fa | an setpoii | nt | | | |
| Po | 00 os1 End | _ | 23456 | [%] 789, | Entry menu "Fan setpoint". A fan speed value is shown. Enter the desired value with the operating button. Setting range: 40 % up to 100 % Press the "Ok" button to confirm. |
| | Close | 0 | k | | |

When entering a value outside the setting range, the message "invalid value" appears. Press the operating button to confirm with "Ok" and repeat the entry with a correct value.

| Fixe | d value | | | | | |
|------------------------|------------------------|---------|---|----------|---------------|--------------------------------|
| \ Te | \ Temperature setpoint | | | | | |
| No. | Value | Day | H:M:S | | Activity | |
| 1 | 35.00 | Monday | 08:00 | 0:00 | Active | |
| 2 | 40.00 | Monday | 10:00 | 0:00 | Active | |
| 3 | 35.00 | Monday | 18:0 | 0:00 | Active | Week program editor (example). |
| 4 | 10.00 | Monday | 20:0 | 0:00 | Active | Press the "Menu" button. |
| 5 | 35.00 | Tuesday | 08:00 | 0:00 | Active | |
| 6 | 40.00 | Tuesday | 10:0 | 0:00 | Active | |
| 7 | 35.00 | Tuesday | 18:0 | 0:00 | Active | |
| 8 | 10.00 | Tuesday | 20:0 | 0:00 | Active | |
| | \bigtriangleup | Mer | u | ~ | $\overline{}$ | |
| | | | | | | 1 |
| Fixe | d value | | 0 | 8.03.201 | 5 05:05:06 | |
| | ogram m | | | | | |
| Sele | ct parame | ter | | | _ | |
| Save | e and exit | | | | | |
| Save | and run | | Submonu "Drogrom mon" | | | |
| Exit (without saving!) | | | | | | Submenu "Program menu". |
| | | | Select the desired function and press the operating button. | | | |

10.3.8 Saving the week program and leaving the program editor

With the "Close" button, The controller returns to the week program editor.

| Select parameter | Changing between temperature and fan speed. |
|------------------------|---|
| Save and exit | Saving the program. The controller returns to the "Week programs" submenu. You can now select and start the week program as described in chap. 10.1. |
| Save and run | Saving the program and start it. Enter the start date and time, see chap. 10.1. If another time or week program is running, the program is only saved, but not started. A corresponding message is displayed. |
| Exit (without saving!) | Attention: the program is not saved. After a security question the con- troller returns to the initial view. |
| Button "Close" | The controller returns to the week program editor. You can continue programming. |

Close

With "Exit (without saving!)" you exit the program editor without saving the program. There is a security question first:

| Fixed value | 08.03.2015 05:05:06 | |
|-------------------------------|---------------------|---|
| \ Exit (without saving) | | |
| Do not exit | | |
| Really exit (without saving!) | | Submenu "Exit (without saving)". This a security question. Select the desired function and press the operating button. Then there is an additional security question. Select the desired function and press the operating button. |
| Close | | |

If you selected "Really exit (without saving!)", the controller goes back to the initial view.

10.4 Deleting a week program

To delete a week program, go to *Menu > Programs > Week program > Delete* or *Menu > Programs > Week program > Delete all*

| Fixed value | 08.03.2015 05:05:06 | |
|-------------------------|---------------------|---------------------------------|
| \ Programs\Week program | | |
| Start | | |
| Stop | | |
| Edit | | Submenu "Week program". |
| Create | | Select "Delete" or "Delete all" |
| Rename | | and press the operating button. |
| Delete | | |
| Delete all | | |
| | | |
| Close | Home | |

If you selected "Delete", select then the week program to be deleted and press the operating button.

If you selected "Delete all", all week programs will be deleted in the controller.

Before deleting there is a security question:

| Fixed value \ Confirm operation Do not delete Really delete | 08.03.2015 05:05:06 | Submenu "Confirm operation". This a security question. Select the desired function and press the operating button. |
|--|---------------------|--|
| Close | Home | |

Go back to the initial view with "Home".

11. Key lock

The key lock function serves to block the access to the controller. When the "key lock" is activated, the controller remains in the actual view and can only be changed when entering the current password.

To configure the key lock function, go to *User > Key lock*

| Fixed value \ User | 08.03.2015 05:05:06 | |
|---------------------------|---------------------|---|
| Key lock | | |
| Show event list | | |
| | | "User" menu. Select "Key lock" and press the operating button. |
| Close | Home | |
| P' 1 - 1 | | |
| Fixed value \ Key lock | 08.03.2015 05:05:06 | |
| Key lock On | | |
| Automatic key lock | | |
| Password | | Submenu "Key lock". Select the desired function and press the operating button. |
| Close | Home | |
| | Tierrie | |

| Key lock On | The key lock is directly activated |
|--------------------|--|
| Automatic key lock | The key lock is activated automatically after a defined waiting time |
| Password | Change password for unlocking. Factory setting: 0000 |

11.1 Directly activating the key lock function

To directly activate the key lock, go to User > Key lock > Key lock On

| Fixed value \ Key lock Key lock On Automatic key lock Password | 08.03.2015 05:05:06 | Submenu "Key lock". Select "Key lock On" to activate the key lock function immediately and press the operating button. |
|--|---------------------|---|
| Close | Home | |

--0

This symbol on the controller display indicates that the "key lock" function is activated.

The controller remains in the initial view and may be operated only after entering the current password.

| Enter password | 08.03.2015 05:05:06 | |
|----------------|---------------------|-------------------|
| \ Key lock pas | sword | |
| | | |
| _ | | |
| | | |
| | | |
| | | |
| | | |
| Ok: A B C D I | ef () | 1 2 3 4 5 6 7 8 9 |
| | | |
| Close | Ok | Home |

Entry menu "Key lock password". Enter the desired password with the operating button. Factory setting is 0000 Press the "Ok" button to confirm.

11.2 Automatic key lock

To configure the automatic key lock, go to User > Key lock > Automatic key lock

| Fixed value | 08.03.2015 | 05:05:06 | |
|----------------------|------------|----------|--|
| \ Automatic key lock | | | |
| Automatic key lock | | | |
| Waiting time [min] | | | |
| | | | Submenu "Automatic Key lock" Select the desired function and press the operating button. |
| Close | Ho | me | |



Under "Waiting time [min]" you can enter the waiting time, after which the key lock will be automatically activated. This time starts running off after the last entry to the controller. To enter it, go to User > Key lock > Automatic key lock > Waiting time [min]

| Enter wait time | [min] 08 | 3.03.2015 05:05:06 | |
|-----------------|----------|--------------------|---|
| \ User\Waiting | time | | |
| Pos1 End Of | < 0 1 23 | 456789 | Entry menu "Waiting time". Enter the desired interval with the operating button. This interval starts running off after the last action on the controller. If the automatic key lock function is enabled, it will become active after this time. Factory setting: 1 minute. Press the "Ok" button to confirm. |
| Close | Ok | Home | |

To activate the automatic key lock function with the pre-configured waiting time, select User > Key lock > Automatic key lock > Automatic key lock

| Fixed value | | 08.03.2015 | 05:05:06 | |
|----------------|----------------|------------|----------|---|
| \ Automatic ke | ey lock On/Off | | | |
| Automatic key | lock : On | | | Submenu "Automatic Key lock On/Off". |
| | | | | The current setting is displayed. |
| | | | | To change the setting, press the operating button. |
| | | | | "Automatic Key lock : On" = Automatic key lock func- tion activated. The time set under "Waiting time" be- gins running off |
| | | | | "Automatic Key lock : Off" = Automatic key lock func- tion deactivated |
| Close | | Но | me | |

Now the waiting time starts running off.

Go back to the initial view with "Home".

| Fixed value | | 0 | 8.03.2015 05:0 | 5:06 |
|------------------------|---------|--------------|----------------|-----------|
| | Setp | oint | Actual val | ue |
| Temperature [Fan [| | 25.0)0.0 | | h |
| | | | | A Io |
| ▶• ◀ - • | | | | |
| User | Quick n | nenu | Menu | |

Initial view.

As soon as the waiting period has expired, the "key lock" symbol is displayed.

The controller remains in the initial view and may be operated only after entering the current password.

After further entries to the controller, the waiting period begins running again, since the automatic keylock function remains active until turning it off manually.



11.3 Changing the password for unlocking the key lock

To change the password for unlocking the key lock, go to User > Key lock > Password


| Fixed value 08.03.2015 | 05:05:06 |
|-----------------------------------|--|
| \ Key lock | |
| Key lock On | |
| Automatic key lock | |
| Password | |
| | Submenu "Key lock". |
| | Select "Password" |
| | and press the operating button. |
| | |
| | |
| | |
| Close Ho | me |
| A accurity question is displayed: | |
| A security question is displayed: | |
| Fixed value 08.03.2015 | 05:05:06 |
| \ Key lock \ Change password | |
| Do not change | |
| Change password | |
| | Submenu "Change password". |
| | To change the password, |
| | select "Change password" |
| | and press the operating button. |
| | |
| | |
| | |
| Close Ho | me |
| | |
| Key lock password 08.03.2015 | 05:05:06 |
| \ Current password | |
| | |
| - | |
| | Entry menu "Current password". |
| | Enter the desired password with the operating but ton. Factory setting is 0000 |
| | Press the "Ok" button to confirm. |
| | |
| Ok: A B C D E F 0 123456 | 789 |
| | |
| | ome |
| Close Ok Ho | |
| | d modification. Without the correct password, unlocking the |
| | a mounication. Without the correct password, unlocking the |

key lock is impossible.

12. General controller settings

In the "Settings" submenu, you can enter the date and time, select the language for the menus and the desired temperature unit, perform the configuration for the controller's communication functions, and reset the controller to factory settings.

To access the "Settings" submenu, go to *Menu* > Settings

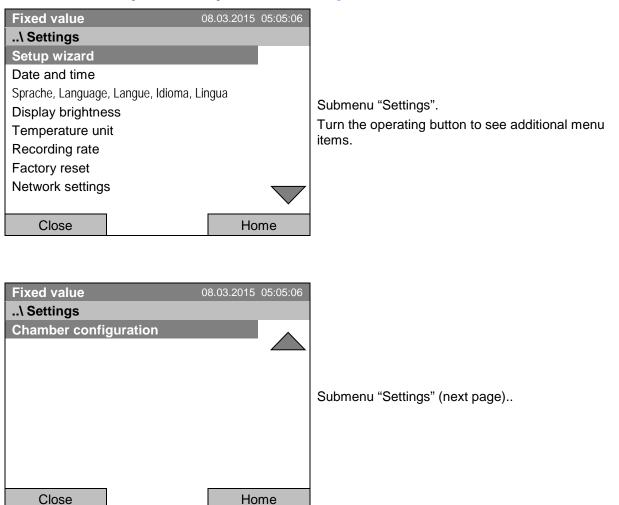


| Setup wizard | Chap. 12.1 |
|--|---|
| Date and time | Setting date and time, chap. 12.2 |
| Sprache, Language, Langue, Idioma, Lingua | Selecting the controller's menu language, chap. 12.3 |
| Display brightness | Adjusting display brightness by turning the operating button |
| Temperature unit | Selecting the temperature unit, chap. 12.5 |
| Recording rate | Defining the recording rate for data storage, chap. 12.6 |
| Factory reset | Factory reset, chap. 12.7 |
| Network settings | Network configuration, chap. 12.8 |
| Chamber configuration | Menu for service purpose – Display and entry of chamber data (serial no., special application no., option object temperature, chap. 12.10 |

12.1 Setup wizard

The setup wizard will guide you sequentially through the important menus to configure your chamber

- Menu language
- Chamber name
- Date and time
- IP address
- Subnet mask
- Network name
- Gateway
- DNS 1
- DNS 2

Then the controller returns to the initial view.

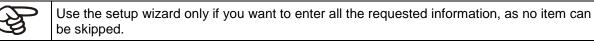


You can configure the network settings (IP address, and the following) only if the DHCP status is Off, otherwise the DHCP server would assign the network configuration.

If you try configuring any network settings while DHCP is enabled, (i.e. settings from menu item "IP Address" on in the setup wizard), the message "DHCP enabled!" is shown. After confirming with Ok the Setup wizard is cancelled and the controller returns to the initial display. The settings made up to then remain effective.

12.2 Date and time settings

To access the date and time settings, go to Menu > Settings > Date and time

| Fixed value | 08.03.201 | 5 05:05:06 | |
|-----------------|-----------|--------------------|---|
| \ Date and time | e | | |
| Set date | | | |
| Set time | | | |
| | | | Submenu "Date and time". |
| | | | Select the desired function and press the operating but |
| | | \bigtriangledown | |
| | | | |
| Close | H | ome | |

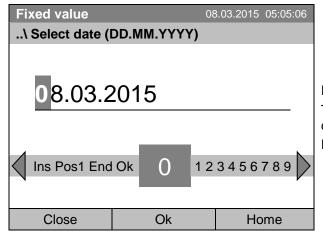

Entry menu "Select date".

The current date is shown. If it is incorrect, enter the correct date with the operating button. Press the "Ok" button to confirm.

Function "Set time"

Function "Set date"

| Fixed value | | 3.03.2015 05:05:06 | |
|--------------|-----------|--------------------|---|
| | | | |
| 05:05:0 |)6 | | Entry menu "Set time". The current time is shown. If it is incorrect, enter the correct time with the operating button. |
| Ins Pos1 End | Ok () 123 | 3 4 5 6 7 8 9 | Press the "Ok" button to confirm. |
| Close | Ok | Home | |

There is no automatic switch to local light-saving times because this could lead to problems with data seeming to be missing or being overwritten in the data base.

Go back to the "Settings" menu with "Close" or to the initial view with "Home".

12.3 Selecting the menu language of the T4.12 controller

The T4.12 chamber controller communicates via a comprehensible menu navigation in plain text in a selectable language.

| wenu > Settings | > Spracne, Language | , Langue, | iaioma, |
|------------------|-----------------------------|-------------|-----------------|
| Fixed value | 08.03.20 | 15 05:05:06 | |
| \ Sprache, Langu | iage, Langue, Idioma, Lingi | ua | |
| Deutsch | | | |
| English | | | |
| Français | | | Quiltant |
| Español | | | Subm |
| Italiano | | | Select and p |
| | | | and p |
| | | | |
| | | | |
| | | | |
| Close | H | Home | |

To select the desired menu language, go to *Menu* > *Settings* > *Sprache, Language, Langue, Idioma, Lingua*

> Submenu "Language". Select the desired language and press the operating button.

If you selected the menu language, the controller returns to the "Settings" menu.. Otherwise, go back to the "Settings" menu with "Close" or to the initial view with "Home".

12.4 Setting display brightness

To select the display brightness, go to Menu > Settings > Display brightness

| Fixed value\ Display brigi | | 8.03.2015 | 05:05:06 | |
|----------------------------|------|-----------|----------|---|
| | | | | Submenu "Display brightness". Select the desired setting by turning the operating button. Confirm by selecting "Save". |
| Close | Save | Но | me | |

Go back to the "Settings" menu with "Close" or confirm the change with "Save": The controller returns to the initial view.

12.5 Changing the temperature unit

To select the temperature unit, go to Menu > Settings > Temperature unit

| Fixed value | 08.0 | 03.2015 | 05:05:06 | |
|-----------------|------|---------|----------|---|
| \ Temperature | unit | | | |
| Celsius [°C] | | | | |
| Fahrenheit [°F] | | | | |
| Kelvin [K] | | | | Submenu "Temperature unit". |
| | | | | Select the desired temperature unit" and press the operating button. |
| | _ | | | |
| Close | | Ho | me | |

If you selected the temperature unit, the controller returns to the "Settings" menu.

Otherwise, go back to the "Settings" menu with "Close" or to the initial view with "Home".

12.6 Defining the data recording rate

To define the recording rate for data storage on the SD card, go to Menu > Settings > Recording rate

| Fixed value | 30 | 3.03.2015 05:05:06 | |
|----------------------------|---------|--------------------|---|
| \ Recording rate (seconds) | | | |
| 60 | 5 () 78 | 9. – Del Ins | Entry menu "Recording rate". The current recording rate is shown. Enter the de- sired value with the operating button. Shortest possible value: 60 seconds. Press the "Ok" button to confirm. |
| Close | Ok | Home | |

When entering a value smaller than 60 seconds, the message "invalid value" appears. Press the operating button to confirm with "Ok" and repeat the entry with a correct value.

Go back to the "Settings" menu with "Close" or to the initial view with "Home".

Note:

After a period of 9 years, the controller starts overwriting the oldest values on the storage medium. This is independent of the selected storage interval and the actual operating time of the chamber. In any case, the data can be read out at any time using the function "Export to USB drive" (chap. 13.1) and stored externally.




12.7 Factory reset

The "factory reset" function allows resetting the controller configuration to the factory settings.

| Risk of data loss! When resetting to factory settings, all controller settings, user data and passwords, will be deleted. Any entered programs, the event list and stored measuring data are NOT affected. |
|---|

To access the "Factory reset" function, go to Menu > Settings > Factory reset

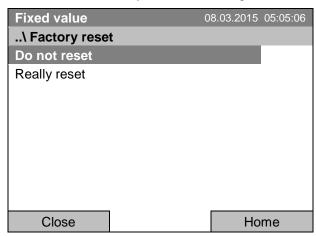

Submenu "Factory reset". This a security question. Select the desired function and press the operating button.

If you selected "Really reset" a message asking to restart the chamber appears. Confirm with "OK". The controller returns to the initial view.

Otherwise, go back to the "Settings" menu with "Close" or to the initial view choosing "Do not reset" or "Home".

12.8 Network configuration

The settings of this submenu are required for networking chambers with an Ethernet interface, e.g. to connect them with BINDER's APT-COM[™] 4 Multi Management Software.

You can display the chamber's IP address that has been assigned by your DHCP server or manually assign the IP address. All necessary configurations for networking the chamber are available in this menu.

To view and configure the network settings, go to Menu > Settings > Network settings

To configure the network settings, the DHCP state must be set to off. Otherwise, the DHCPserver would assign the network configuration.

If you try to configure the network settings with enabled DHCP state, the message "DHCP enabled!" is shown. Confirm with "Ok" to return to the "Network settings" menu.

| Fixed value | 08.03.2015 | 05:05:06 | |
|-----------------------|------------|--------------------|---|
| \ Network settings | | | |
| Show network settings | | | |
| DHCP On/Off | | - | |
| MAC address | | | Output and "National a stitute" |
| IP address | | | Submenu "Network settings". |
| Subnet mask | | | Turn the operating button to see additional menu items. |
| Chamber name | | | items. |
| Network name | | | |
| Gateway | | \bigtriangledown | |
| Close | Но | me | |

BINDER

| Fixed value | 30 | 8.03.2015 05:05:06 | |
|----------------|------|--------------------|---|
| \ Network sett | ings | | |
| DNS 1 | | | |
| DNS 2 | | | • |
| | | | |
| | | | Submany "Notwork acttings" (next nega) |
| | | | Submenu "Network settings" (next page). |
| | | | |
| | | | |
| | | | |
| | - | | |
| Close | | Home | |

| Show network settings | Overview of the entire network configuration |
|-----------------------|--|
| DHCP on/off | Switching on and off the DHCP state |
| MAC address | Display of the chamber's MAC address |
| IP address | Entering the desired IP address |
| Subnet mask | Entering the subnet mask number |
| Chamber name | Entering the name of the cooling incubator |
| Network name | Entering the network name |
| Gateway | Entering the gateway number |
| DNS 1 | Entering the DNS 1 number |
| DNS 2 | Entering the DNS 2 number |

Set the DHCP State on/off:


| Fixed value | 22.08.2012 | 05:05:06 | |
|---------------|------------|----------|---|
| \ DHCP On/Off | | | |
| DHCP : On | | | |
| | | | Submenu "DHCP On/Off". |
| | | | The current DHCP state is displayed. Press the operating button to change it. |
| | | | Then the new DHCP state is displayed |
| | | | "DHCP : On" = DHCP state activated "DHCP : Off" = DHCP state deactivated |
| | | | |
| Close | Но | me | |

Go back to the "Network settings" menu with "Close" or to the initial view with "Home".

You can only configure the following network settings if the DHCP state is set to "Off".

Display the MAC address

To identify the chamber in the Ethernet network you can display the chamber's MAC address.

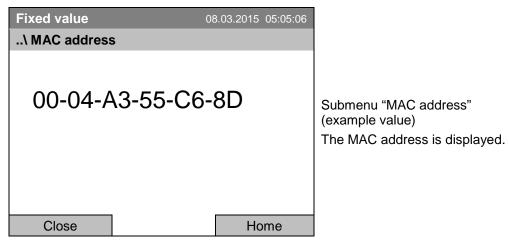

Go back to the "Network settings" menu with "Close" or to the initial view with "Home".

Enter the IP address:

| Fixed value | 22 | 2.08.2012 0 | 5:05:06 | |
|-------------------------|----|--|---------|--|
| \ IP address (n.n.n.n.) | | | | |
| 1 92.168.0.100 | | Entry menu "IP address" (example value) Enter the desired IP address with the operating but ton. Press the "Ok" button to confirm. | | |
| Close | Ok | Hom | ie | |

Go back to the "Network settings" menu with "Close" or to the initial view with "Home".

Enter the subnet mask:

| Fixed value | 22 | 2.08.2012 05:05:06 | | |
|--|----|--------------------|--|--|
| \ Subnet mask (n.n.n.n.) | | | | |
| .\ Subnet mask (n.n.n.) 2 55.255.255.0 Ins Pos1 End Ok 0 1 2 3456789,- | | 56789 - | Entry menu "Subnet mask" (example value) Enter the desired net mask with the operating but- ton. Press the "Ok" button to confirm. | |
| N | | | | |
| Close | Ok | Home | | |

Go back to the "Network settings" menu with "Close" or to the initial view with "Home".

Enter the chamber name:

| Fixed value | 22 | 2.08.2012 05:05:06 | |
|---------------|----|--------------------|---|
| \ Chamber nar | ne | | |
| KB_E6 | | INOPQR | Entry menu "Chamber name" (example) Enter the desired chamber name with the operating button. Press the "Ok" button to confirm. |
| Close | Ok | Home | |

Go back to the "Network settings" menu with "Close" or to the initial view with "Home".

Enter the network name:

| Fixed value | 22 | 2.08.2012 05:05:06 | |
|---------------|----------|--------------------|---|
| \ Network nam | ne | | |
| | 、 | | Entry menu "Network name" |
| K B_E6 | | | (example) |
| | | | Enter the desired network name with the operating button. |
| CDEFGH | IJ K LN | INOPQR | Press the "Ok" button to confirm. |
| Close | Ok | Home | |

Go back to the "Network settings" menu with "Close" or to the initial view with "Home".

Enter the default gateway:

| Fixed value | 22 | 2.08.2012 05:05:06 | |
|----------------|----------|--------------------|---|
| \ Gateway (n.r | n.n.n.) | | |
| 192.168.0.1 | | | Entry menu "Gateway" (example value) Enter the desired gateway with the operating button. |
| Pos1 End O | k 0 1 23 | 456789, | Press the "Ok" button to confirm. |
| Close | Ok | Home | |

Go back to the "Network settings" menu with "Close" or to the initial view with "Home".

Enter the DNS 1 or DNS 2:

| Fixed value | 2 | 2.08.2012 05:05:00 | |
|-----------------|------|--------------------|--|
| \ DNS 1 (n.n.n. | .n.) | | |
| 192.16 | _ | 456789, | Entry menu "DNS 1" or "DNS 2" (example value) Enter the desired number with the operating button. Press the "Ok" button to confirm. |
| Close | Ok | Home | |

Go back to the "Network settings" menu with "Close" or to the initial view with "Home".

12.9 Display of the entire network configuration

To access the overview of the complete network configuration, go to *Menu* > *Settings* > *Network settings* > *Show network settings*

| Fixed value | 08.03.2015 05:05:06 | | |
|---------------|---------------------|---------------|---|
| \ Show networ | k settings | | |
| DHCP | Off | | |
| MAC address | 00-04 | 4-A3-55-C6-8D | |
| IP address | 192.1 | 168.0.100 | |
| Net mask | 255.2 | 255.255.0 | Overview of the network configuration (sample values) |
| Gateway | 192.1 | 168.0.1 | |
| DNS1 | 192.1 | 168.0.1 | |
| DNS2 | 0.0.0 | .0 | |
| Chamber name | KB_E | Ξ6 | |
| BIOS name | KB_E6 | | |
| Close | | Home | |

Go back to the "Network settings" menu with "Close" or to the initial view with "Home".

12.10 Display and entry of the chamber configuration – for service purpose

Information about the chamber, such as chamber type, name, serial number, firmware version, etc. can be viewed under *Menu* > *System information* (chap. 6.4).

To access the device configuration menu, go to Menu > Settings > Chamber configuration

This menu is password protected and only intended for service purposes.

13. Data transfer via USB interface

The USB port is located in the instrument box.

To access the submenus for data transfer, go to Menu > Import/Export

| Fixed value | 08.03.2015 05:05:06 | |
|------------------------------|---------------------|---|
| \ Import/Export to USB drive | e | |
| Export to USB drive | | |
| Import from USB drive | | |
| | | Submenu "Import/Export to USB drive". |
| | | Select the desired type of communication via USB port (data export or import) and press the operating button. |
| Close | Home | |
| It is possible that sor | ne USB devices are | not recognized due to compatibility issues. In this |

Exporting data to USB drive


S

13.1

Insert the USB stick or the plug of your USB drive into the USB port in the instrument box.

case, please use a USB storage device from a different manufacturer.

To configure data export to USB media, go to Menu > Import/Export > Export to USB drive

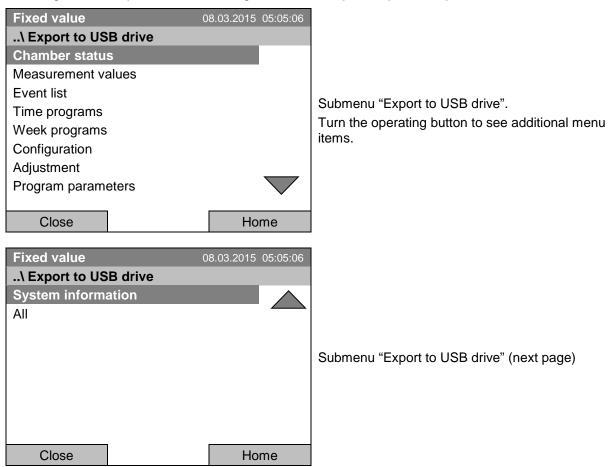



| Chamber status | Actual chamber status, including operating mode, set-points etc. | |
|--------------------|--|--|
| Measurement values | Measured data | |
| Event list | List of status information and errors (see chap. 15) | |
| Time programs | All stored time programs | |
| Week programs | All stored week programs | |
| Configuration | (Service only) | |
| Adjustment | Adjustment data | |
| Program parameters | (Service only) | |
| System information | (Service only) | |
| All | All data | |

Select the desired data type and press the operating button. Data will be written to the connected media.



If no USB device has been connected, the message "No USB device found" is displayed. It disappears after inserting the USB stick or the plug of your USB drive into the USB port in the instrument box.




This symbol on the controller display indicates that data are being transmitted via the USB port.

13.2 Importing data from USB drive

Insert the USB stick or the plug of your USB drive into the USB port in the instrument box.

To configure data import from USB media, go to *Menu > Import/Export > Import from USB drive*

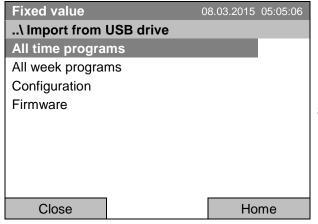

Submenu "Import from USB drive".

Select the desired data type and press the operating button. Data will be imported from the connected media.

| All time programs | All stored time programs | |
|-------------------|--------------------------|--|
| All week programs | All stored week programs | |
| Configuration | (Service only) | |
| Firmware | (Service only) | |



If no USB device has been connected, the message "No USB device found" is displayed. It disappears after inserting the USB stick or the plug of your USB drive into the USB port in the instrument box.



This symbol on the controller display indicates that data are being transmitted via the USB port.

14. Notifications and Alarms

14.1 Notifications overview

| lcon | Signification | | lcon | Signification |
|------|------------------------------------|-----------------|---|--|
| ▶•◀ | Fixed value operation | | <u> </u> | Heating active |
| Ċ | Time program operation | g modes | <u>22</u> | Door heating active (KB / KB-UL 240 / 400 / 720) |
| \$ | Time program interrupted | Operating modes | * | Refrigeration active |
| ß | Week program operation | | 0 | Key lock activated |
| Ð | Interior socket (option) turned on | | USB | Copying data via USB |
| | | | $\begin{vmatrix} 1 \bullet \\ 2 \bullet \\ 3 \end{vmatrix}$ | Switching state of the optional zero- voltage relay control outputs (example: control outputs 1 + 2 switched on) |

14.2 Alarms overview

| Icon | Alarm message | Signification | |
|------------|---------------------------------------|---|--|
| ∞ | Safety controller overtemperature | Overtemperature safety controller alarm (class 3.1): selected value of the safety controller exceeded | |
| ₽Ţ! | Safety controller overtemperature | Overtemperature safety controller alarm with option class 3.3: selected value of the safety controller exceeded | |
| J. | Safety controller undertemperature | Undertemperature safety controller alarm with option class 3.3: fallen be- low selected value of the safety controller | |
| ₽ | Temp. range | Tolerance range alarm: Having reached the set-point, the temperature deviates by more than $+/- 2$ °C from the set-point for more than 10 minutes, or the temperature doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door | |



You can activate / deactivate the buzzer in the "Alarms" submenu (chap. 14.5.3).

With an activated buzzer there is an **audible alert** with an alarm. You can reset it in the "Alarms" submenu for alarm acknowledgement pressing the "Reset" button (chap. 14.4). The alarm symbol will only disappear when the cause of the alarm has been remedied.

An optional **zero-voltage relay alarm contact for temperature** (chap. 19.3) is available. It is activated in case of the alarm messages with temperature deviations) as well as in case of a power failure and when turning off the main power switch.

For appropriate actions in the event of an alarm, please refer to chap. 21.3 "Troubleshooting".

14.3 Alarm status

An alarm message can appear in 3 different states:

Status "set"


- Active alarm.
- The corresponding alarm icon is displayed in the initial view The buzzer sounds (if activated).
- The "Info" button in the initial view leads to the "Alarms" submenu for alarm acknowledgement.
- Press the "Reset" button in the "Alarms" submenu for alarm acknowledgement to mute the buzzer and confirm the alarm.

Status "acknowledged"

- Active alarm.
- The alarm was acknowledged.. The buzzer is off.
- The alarm cause is still valid. Therefore, the alarm icon remains displayed in the initial view.
- The alarm message figures in the list of active alarms.

Status "cleared"

- The alarm cause has disappeared.
- · The alarm icon is not displayed any longer
- The alarm message has disappeared from the list of active alarms.
- The alarm message remains in the Event list for information.

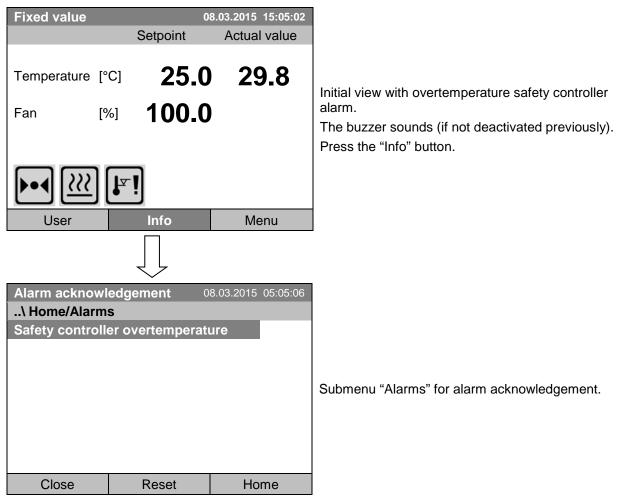

14.4 Confirming a "set" alarm

Press the "Reset" button to confirm the alarm. If there is another active "set" alarm, you can also confirm it with the "Reset" button. After confirming all active "set" alarms, the buzzer is off, and the initial view is displayed.

As long as an alarm is still active, i.e. as long as the alarm cause is still valid, the alarm icon continues to be displayed in the initial view.

14.5 Alarm configuration and overview

To access the alarm lists and configuration menu, go to Menu > Alarms

| Fixed value | 0{ | 8.03.2015 | 05:05:06 | |
|--|----|-----------|----------|--|
| Active alarms | | | | |
| History Buzzer test Buzzer On/Off Alarms On/Off | | | | Submenu "Alarm". Select the desired function and press the operating button. |
| Close | | Но | me | |

| Active Alarms | List of the active alarms (status "set" or "acknowledged"). | | |
|--|--|--|--|
| History | List of all alarms (status "set" or "acknowledged" or "cleared"). | | |
| Buzzer test | Testing the alarm buzzer, chap. 14.5.3 | | |
| Buzzer On/Off Activating / deactivating the alarm buzzer, chap. 14.5.3 | | | |
| Alarms On/Off | Activating / deactivating the alarm functions. Off: Alarm buzzer off, no alarm icons displayed. No alarm messages will be displayed. | | |

14.5.1 List of active alarms

To access the overview list of active alarms, go to Menu > Alarms > Active alarms

| Fixed value | 0 | 8.03.2015 | 05:05:06 | |
|------------------|------------------|-----------|----------|--|
| \ Active alarm | s | | | |
| Safety controlle | er overtemperati | ure | | |
| | | | | |
| | | | | Submenu "Active alarms". |
| | | | | All active alarms with status "set" or "acknowledged", |
| | | | | are listed. |
| | | | | |
| | | | | |
| | 1 | | | |
| Close | | Но | me | |

If no alarm is active ("set" or "acknowledged"), no message will be displayed in this window.

14.5.2 History - list of all alarms

To access the overview list of all alarms, go to Menu > Alarms > History

This list indicates the moment when an alarm was set and when cleared.

| Fixed value 08. | 03.2015 15:05:06 | |
|---|------------------|--|
| \ History | | |
| 08.03.2015 : Messages of indicate | ed day | |
| 14:39:48 Alarm set Safety controller overte | emperature | Submenu "History" (example). |
| 15:03:22 Alarm cleared Safety controller ov | vertemperature | The list shows when the alarms of the current day were triggered and when cleared. The most recent message appears at the end of the list. |
| | | The information that an alarm has been acknowl- edged is shown in the event list. |
| Close | Home | |

When there is more information than one page, you can scroll the list in both directions with the operating button.

To select a different date, select "Messages of indicated day" and press the operating button. You can enter the desired date through an entry menu.

| Fixed value | | 08.03.2015 | 05:05:06 |
|------------------|-------------|------------|---|
| \ Select date (I | DD.MM.YYYY) | | |
| | | | |
| 08.03.2 | 2015 | | |
| | | | |
| 1 | | | |
| Ins Pos1 End | Ok () ' | 123456 | 789 🔪 |
| | | | r |
| Close | Ok | Ho | ome |
| 0.000 | ÖK | 110 | ,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,, |

Entry menu "Select date". The current date is shown. Enter the desired date with the operating button. Press the "Ok" button to confirm.

The alarm list of the selected date is displayed.

The entire sequence of the alarm events (set – acknowledged – cleared) is shown in the event list (chap. 15).

To access the event list, go to *Menu > Event list* or *User > View event list*

| Fixed value | 08.03.2015 15:05:06 | |
|---------------------------------------|--------------------------------|--|
| \ Event list | | |
| 08.03.2015 : Messages of | f indicated day | |
| 14:17:20 Fan setpoint 50.00 |) | |
| 14:35:12 Fan setpoint 100.0 | 00 | Submenu "Event list" (example). |
| 14:39:48 Alarm set Safety contr | oller overtemperature | The events and alarm messages of the current day |
| 14:40:19 Alarm acknowledged S ture | Safety controller overtempera- | are displayed. The most recent message appears at the end of the list. |
| 15:03:22 Alarm cleared Safety of | controller overtemperature | |
| | | |
| | | |
| Close | Home | |

BINDER

14.5.3 Activating, deactivating, and testing the alarm buzzer

Alarm buzzer test

To access the functional test of the alarm buzzer, go to Menu > Alarms > Buzzer test

| Fixed value | 30 | 8.03.2015 05:05:06 | |
|-----------------|----|--------------------|--|
| \ Buzzer test | | | |
| Buzzer test: Of | f | | Submenu "Buzzer test". |
| | | | The current setting is displayed. |
| | | | Press the operating button to turn on or off the buzzer for test purpose. |
| | | | The modified setting is displayed |
| | | | "Buzzer test: On" = Buzzer turns on "Buzzer test: Off" = Buzzer turns off |
| Close | | Home | |

When turned on, the alarm buzzer emits an intermittent signal. To turn this off, change the test function to "Buzzer test: Off".

Activating / deactivating the alarm buzzer

To activate or deactivate the alarm buzzer, go to Menu > Alarms > Buzzer On/Off

| Fixed value | C | 8.03.2015 | 05:05:06 | |
|-----------------|--------|-----------|----------|--|
| \ Buzzer On/O | ff | | | |
| Buzzer activati | on: On | | | Submenu "Buzzer On/Off". |
| | | | I | The current setting is displayed. |
| | | | | To change the setting, press the operating button. |
| | | | | The modified setting is displayed |
| | | | | |
| | | | | "Buzzer activation: On" = Buzzer will turn on in the event of an alarm |
| | | | | "Buzzer activation: Off" = Buzzer is deactivated |
| | | | | |
| Close | | Но | me | |

14.5.4 Activating / deactivating all alarm functions

To access the alarm settings, go to Menu > Alarms > Alarms On/Off

| Fixed value | 08.03.2015 05:05:06 | |
|--|---------------------|---|
| \ Alarm functions (Alarm functions: Or | | Submenu "Alarm functions On/Off". The current setting is displayed. To change the setting, press the operating button. The modified setting is displayed "Alarm functions: On" = The alarm functions of the controller are active "Alarm functions: Off" = All alarm functions are deac- tivated |
| Close | Home | |

Go back to the initial view with "Home".

15. Event list

The "Event list" displays status information and errors of the current day. You can also access the events of past days.

To access the event list, go to *Menu > Event list* or *User > Show event list*

| 10:35:12 Fan setpoint 80.00 | Fixed valu | e | 08.03.2015 | 15:05:06 | |
|--|------------------------------------|---------------------|-------------|----------|--|
| 10:11:59Power supply was interrupted.Submenu "Event list" (example).10:12:05Common controller mode OFFThe events of the current day are displayed. The most recent message appears at the end of the list10:35:12Fan setpoint 80.00To acknowledge an activated alarm, proceed as of | \ Event lis | st | | | |
| 10:12:05Common controller mode OFFThe events of the current day are displayed. The most recent message appears at the end of the lis10:17:20Common controller mode ONmost recent message appears at the end of the lis10:35:12Fan setpoint 80.00To acknowledge an activated alarm, proceed as determined and the list | 10:11:49 | Time program sto | oped | | |
| 10:17:20Common controller mode ONmost recent message appears at the end of the list10:35:12Fan setpoint 80.00To acknowledge an activated alarm, proceed as d | 10:11:59 | Power supply was i | nterrupted. | | Submenu "Event list" (example). |
| 10:35:12Fan setpoint 80.0011:04:12Fan setpoint 100.00To acknowledge an activated alarm, proceed as d | 10:12:05 | Common controller | mode OFF | - | |
| 11:04:12 Fan setpoint 100.00 To acknowledge an activated alarm, proceed as d | 10:17:20 Common controller mode ON | | | | most recent message appears at the end of the list. |
| | 10:35:12 | Fan setpoint 80.00 | | | |
| | 11:04:12 | Fan setpoint 100.00 |) | | To acknowledge an activated alarm, proceed as described in chap. 14.4. |
| | | | | | |
| Close Reset Home | Close | Reset | Ho | me | |

Turn the operating button to see additional data.

| Fixed valu | | | | | | | | |
|------------|------|---------------------------|----------|----|--|--|--|--|
| \ Event li | | | | | | | | |
| 08.03.2015 | Mes | Messages of indicated day | | | | | | |
| 09:01:59 | Usei | r management o | lisabled | | | | | |
| 09:12:05 | Lang | guage loaded | | | | | | |
| 09:17:20 | Fan | setpoint 80.00 | Submenu | | | | | |
| 09:35:12 | Fan | The event | | | | | | |
| 10:04:12 | New | | | | | | | |
| 10:06:46 | Time | | | | | | | |
| | | | | | | | | |
| | | | | · | | | | |
| Close | | Reset | Ho | me | | | | |

Submenu "Event list" (example). The events of the current day are displayed.

When there is more information than one page, you can scroll the event list in both directions with the operating button.

To select a different date, select "Messages of indicated day" and press the operating button. You can enter the desired date through an entry menu.

| Fixed value | 30 | 8.03.2015 05:05:06 | |
|-----------------|-------------|--------------------|---|
| \ Select date (| DD.MM.YYYY) | | |
| 08.03.2 | _ | 3456789 | Entry menu "Select date". The current date is shown. Enter the desired date with the operating button. Press the "Ok" button to confirm. |
| Close | Ok | Home | |

The event list of the selected date is displayed.

16. Graphical display of the measured values

To access the graphical display, go to Menu > Measurement chart

| F | ixe | d value | | | | 08.03 | 3.2015 | 05:05:06 | |
|-------------------|-------------------|---------------------|-----|----|--------|-------|--------|-------------|--|
| | | 0:0 | 0:5 | (| 0:11 | 0:16 | 0 | :22 | |
| Temperature | Objecttemperature | | | | | | | | Measurement chart view. (sample view with optional chamber equipmen Press the "Settings" button. |
| Ou. Ou. Ou. | 2 | Close | | Se | ttings | | Hc | ome | |
| | _ | d value ngs | | | - | 08.03 | 3.2015 | 05:05:06 | |
| S | am | pling ra ay rang | | | | | | | |
| | - | meters | | | | | | | Submenu "Settings". Select the desired function and press the operating button. |
| | | Close | | | | | Ho | me | |


16.1 Setting the sampling rate

To define the sampling rate, go to *Menu > Measurement chart > Settings > Sampling rate*

| Fixed value | 08.03.2015 | 05:05:06 | |
|---------------|------------|----------|---|
| Sampling rate | | _ | |
| Five seconds | | | |
| Ten seconds | | | |
| One minute | | | 0 |
| Five minutes | | | Submenu "Sampling rate". |
| Ten minutes | | | Select the desired interval and press the operating button. |
| | | | and press the operating button. |
| | | | |
| | | | |
| | | | |
| Close | Но | me | |

16.2 Defining the display range

To define the sampling rate, go to *Menu > Measurement chart > Settings > Display range*

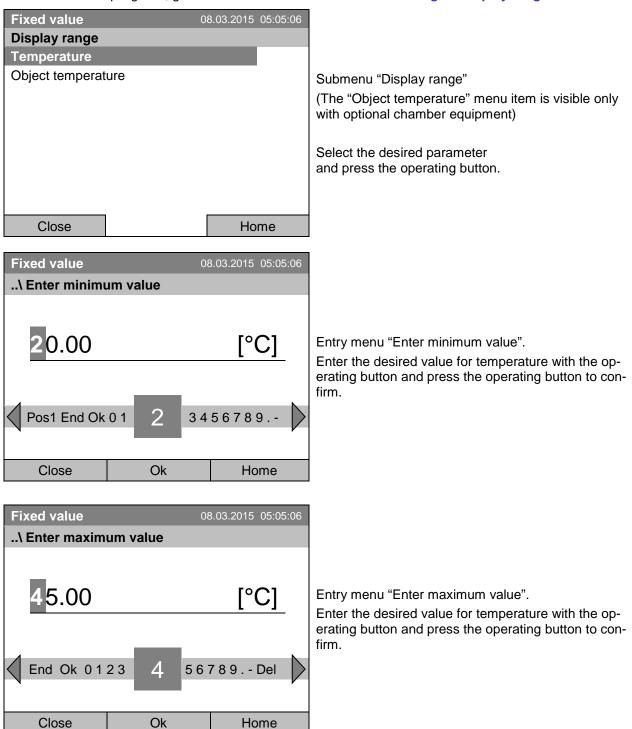

You can now define the view range of other parameters or press the "Close" button twice to go back to the graphic display.

16.3 Selecting the parameters

Here you can select the parameters, which shall be displayed graphically. To select the parameters, go to *Menu > Measurement chart > Settings > Parameters*

| Fixed value | 0 | 8.03.2015 05:05:06 | |
|-------------------|--------|--------------------|---|
| Parameters | | | Submenu "Parameters" |
| Temperature Or | n | | (The menu items "Object temperature" and "Control |
| Object temperatu | ure On | | outputs" are visible only with optional chamber |
| Control outputs (| On | | equipment) |
| | | | |
| | | | The current setting for each parameter is displayed. |
| | | | Select the desired parameter |
| | | | and press the operating button to change its status On/Off. |
| | | | The modified setting is displayed. |
| Close | | Home |] |

Press the button "Close" twice to go back to the graphic display. If any parameter was set to "Off", it will not be included in the graphical display.

17. Temperature safety devices

17.1 Over temperature protective device (class 1)

The chamber is equipped with an internal temperature safety device class 1 acc. to DIN 12880:2007. It serves to protect the chamber and prevents dangerous conditions caused by major defects.

If a temperature of about 110 °C / 230 °F is reached, the over temperature protective device permanently turns off the chamber. The user cannot restart the device again. The protective cut-off device is located internally. Only a service specialist can replace it. Therefore, please contact an authorized service provider or BINDER Service.

17.2 Overtemperature safety controller (temperature safety device class 3.1)

The chamber is regularly equipped with an electronic overtemperature safety controller (temperature safety device class 3.1 according to DIN 12880:2007).

The overtemperature safety controller serves to protect the chamber, its environment and the contents from exceeding the maximum permissible temperature. In the case of an error, it limits the temperature inside the chamber to the entered safety controller set-point.

Please observe the DGUV guidelines 213-850 on safe working in laboratories, issued by the employers' liability insurance association (for Germany).

Set the safety controller set point by approx. 2 °C up to 5 °C above the desired temperature set point. Recommended setting: Set-point type "Offset" with safety controller set-point 2 °C (factory setting).

The safety controller is functionally and electrically independent of the temperature control system. If an error occurs, it performs a regulatory function.



Check the setting regularly and adjust it following changes of the set-point or charge.

The safety controller only activates after the set-point has been reached once.

17.2.1 Safety controller modes

You can select between "Limit (absolute)" and "Offset (relative)" safety controller mode

• Limit: Absolute maximum permitted temperature value

This setting offers high safety as a defined temperature limit will not be exceeded. It is important to adapt the safety controller set-point after each modification of the temperature set-point. Otherwise, the limit could be too high to ensure efficient protection, or, in the opposite case, it could prevent the controller from reaching an entered set-point outside the limit range.

• Offset: Maximum overtemperature above any active temperature set point. The maximum temperature changes internally and automatically with every set-point change.

This setting is recommended for program operation. It is important to check the safety controller setpoint and safety controller mode occasionally, as it does not offer a fix, independent limit temperature value, which would never be exceeded.

Example: Desired temperature value: 40 °C, desired safety controller value: 45 °C.

Possible settings:

| Temperature set point | Safety controller mode | Safety controller set-point |
|-----------------------|------------------------|-----------------------------|
| 40 °C | Limit (absolute) | 45 °C |
| 40 C | Offset (relative) | 5 °C |

17.2.2 Setting the safety controller

To display and to change the current safety controller settings in the "safety controller" submenu, go to *Menu* > *Safety controller*


Safety controller mode: selection between Limit (absolute) and Offset (relative)

| Fixed value | 08 | 8.03.2015 | 05:05:06 | |
|-----------------|-------|-----------|----------|---|
| \ Safety contro | oller | | | |
| Mode | | | | |
| Setpoint | | | | |
| Show settings | | | | Submenu "Safety controller" (view with standard chamber equipment) |
| | | | | Select "Mode" and press the operating button. |
| Close | | Но | me | |

Or



| Fixed value | 08.03.2015 05:05:06 | |
|---------------------|---------------------|--|
| \ Safety controller | | |
| Mode | | |
| Overtemperature | | |
| Undertemperature | | Submenu "Safety controller" |
| Show settings | | (view with optional chamber equipment) |
| | | Select "Mode" |
| | | and press the operating button. |
| | | |
| | | |
| | | |
| Close | Home | |
| | | |

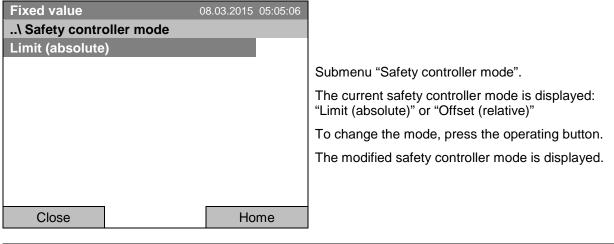

When changing the safety controller mode, the safety controller setpoint which had been active before in this mode becomes active again.

Entering the safety controller setpoint


| Fixed value | 08.03.2015 05:05:06 | |
|---------------------|---------------------|---|
| \ Safety controller | | |
| Mode | | |
| Setpoint | | |
| Show settings | | Submenu "Safety controller" (view with standard chamber equipment) Select "Setpoint" and press the operating button. |
| Close | Home | |

Or

| Fixed value | 08.03.2015 | 05:05:06 | |
|-----------------------------------|------------|----------|--|
| \ Safety controller | | | |
| Mode | | _ | |
| Overtemperature | | | |
| Undertemperature Show settings | | | Submenu "Safety controller" (view with optional chamber equipment) Select "Overtemperature" and press the operating button. |
| Close | Hc | ome | |

In case of a temperature safety device class 3.1 combined with optional chamber equipment, the safety controller setpoint is indicated as "Overtemperature". In this case an additional parameter "Undertemperature" is shown. This latter has no function; if you select it, the message "Setpoints not available" appears.

You can also access this submenu to directly enter the safety controller setpoint via *Quick menu > safety controller setpoint*

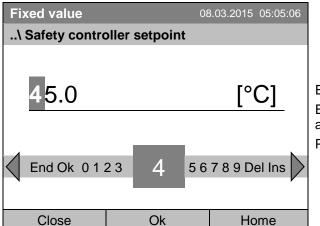

Entry menu "Safety controller setpoint". Enter the desired value with the operating button and press the operating button to confirm. Press the "Ok" button to confirm.

Go one level back with the "Close" button or back to the initial view with "Home".

Overview of the current settings

You can check the current settings of the safety controller:

| Fixed value | 0 | 8.03.2015 | 05:05:06 | |
|-----------------|------|-----------|----------|--|
| \ Safety contro | ller | | | |
| Mode | | | | |
| Setpoints | | | | |
| Show settings | | | | Submenu "Safety controller" (view with standard chamber equipment) Select "Show settings" and press the operating button. |
| Close | | Но | me | |

The overview display shows the set-points and actual values of the main temperature controller and the safety controller and indicates the safety controller mode.

| Fixed value | | 08.03.201 | 5 05:05:06 |
|------------------------|------|-------------|------------|
| | | Setpoint | Actual |
| Chamber temperature | [°C] | 37.0 | 37.6 |
| Safety controller mode | | Limit (abs | olute) |
| Safety controller | [°C] | 40.0 | 37.6 |
| Close | | ŀ | lome |
| Fixed value | | 08.03.201 | 5 05:05:06 |
| | | Setpoint | Actual |
| Chamber temperature | [°C] | 37.0 | 37.6 |
| Safety controller mode | | Offset (rel | ative) |
| Safety controller | [°C] | 3.0 | 37.6 |
| | | | |
| | | | |

Go back to the initial view with "Home".

17.3 Over- and undertemperature safety controller (temperature safety device class 3.3) (option)

With this option the chamber is equipped with an electronic over- and undertemperature safety controller. The combination of overtemperature (class 3.1) and undertemperature (class 3.2) protection is regarded as a safety device class 3.3 acc. to DIN 12880:2007.

This over- and undertemperature safety controller serves to protect the chamber, its environment and the contents from exceeding the maximum or minimum permissible temperature. Please observe the DGUV guidelines 213-850 on safe working in laboratories, issued by the employers' liability insurance association (for Germany). The safety controller is functionally and electrically independent of the temperature control system. If an error occurs, it performs a regulatory function.

The **overtemperature safety controller set-point** is a maximum temperature value, which the chamber will not exceed due to the regulatory function of the safety device. This protection against excessively high temperatures serves to protect the chamber, its environment and the contents. In the case of an error, the safety controller limits the temperature inside the chamber to the entered safety controller set-point.

The **undertemperature safety controller set-point** is a minimum temperature value, which the chamber will not fall below due to the regulatory function of the safety device. This protection against excessively low temperatures serves to protect sensitive loads against under cooling. In the case of an error, the safety controller limits the temperature inside the chamber to the entered safety controller set-point.



Check the setting regularly and adjust it following changes of the set-point or charge.

The safety controller only activates after the set-point has been reached once.

17.3.1 Safety controller modes

You can select between "Limit (absolute)" and "Offset (relative)" safety controller mode

• Limit: Absolute maximum or minimum permitted temperature value

This setting offers high safety as a defined temperature limit will not be exceeded or fallen below. It is important to adapt the safety controller set-point after each modification of the temperature set-point. Otherwise, the limit could be too big to ensure efficient protection, or, in the opposite case, it could prevent the controller from reaching an entered set-point outside the limit range.

• Offset: Maximum permitted over- or under temperature above/ below the active temperature set point. The maximum temperature changes internally and automatically with every set-point change.

This setting is recommended for program operation. It is important to check the safety controller setpoint and safety controller mode occasionally, as it does not offer a fix limit value, which would not be exceeded or fallen below.

Example: Desired temperature value: 40 °C, desired overtemperature safety controller value: 45 °C, desired undertemperature safety controller value: 30 °C.

Temperature
set pointSafety controller
modeOvertemperature safety
controller set-pointUndertemperature safety
controller set-point40 °CLimit (absolute)45 °C30 °C40 °COffset (relative)5 °C10 °C

Possible settings:

17.3.2 Setting the Over- and undertemperature safety controller

To display and to change the current safety controller settings in the "safety controller" submenu, go to *Menu* > *Safety controller*

| Fixed value | 08.03.2015 | 05:05:06 | |
|--|------------|----------|---|
| \ Safety controller | | | |
| Mode | | | |
| Overtemperature Undertemperature Show settings | | | Submenu "Safety controller". Select "Mode" and press the operating button. |
| Close | Ho | ome | |
| Fixed value | 08.03.2015 | 05:05:06 | |
| \ Safety controller mode | | | |
| Limit (absolute) | | | |
| | | | Submenu "Safety controller mode". |
| | | | The current safety controller mode is displayed: "Limit (absolute)" or "Offset (relative)" |
| | | | To change the mode, press the operating button. |
| | | | The modified safety controller mode is displayed. |
| | | | |
| Close | Ho | ome | |

Selection between Limit (absolute) and Offset (relative) safety controller mode



When changing the safety controller mode, the safety controller setpoint which had been active before in this mode becomes active again.

Entering the safety controller setpoint for overtemperature

| Fixed value | 0 | 8.03.2015 05:05:06 | |
|----------------------------------|------|--------------------|--|
| \ Safety contro | ller | | |
| Mode | | | |
| Overtemperatur | е | | |
| Undertemperatur Show settings | e | | Submenu "Safety controller" Select "Overtemperature" and press the operating button. |
| Close | | Home | 1 |

| Fixed value | 30 | 3.03.2015 05:05:06 | |
|----------------|--------------|--------------------|---|
| \ Setpoint ove | rtemperature | | |
| 45.0 | _ | [°C] | Entry menu "Setpoint overtemperature". Enter the desired value with the operating button. Press the "Ok" button to confirm. |
| End , - 012 | 23 4 56 | 7 8 9 Ins Del | |
| Close | Ok | Home | |

You can also access this submenu to directly enter the overtemperature safety controller setpoint via *Quick menu > Safety controller setpoint*

| Fixed value | 08.03.2015 05:05:06 | |
|-----------------------------|---------------------|--|
| \ Safety controller | | |
| Mode | | |
| Overtemperature | | |
| Undertemperature | | Submenu "Safety controller" |
| Show settings | | • |
| | | Select "Undertemperature" with the operating button and press the operating button to confirm. |
| | | |
| | | |
| | | |
| Close | Home | |
| Fixed value | 08.03.2015 05:05:06 | |
| \ Setpoint undertemperature | | |
| | - | |
| - | | |
| 20.0 | [°C] | Entry menu "Setpoint undertemperature". |
| | <u> </u> | Enter the desired value with the operating button. |
| | | Press the "Ok" button to confirm. |
| Pos1 End , - 0 1 2 3 | 456789 Ins | |
| | | |
| Close Ok | Home | |
| Close Ok | поше | |

Entering the safety controller setpoint for undertemperature

Go one level back with the "Close" button or back to the initial view with "Home".

Overview of the current settings

You can check the current settings of the safety controller:

| Fixed value | 08.03.2015 05:05:06 |
|---------------------|---------------------|
| \ Safety controller | |
| Mode | |
| Over temperature | |
| Under temperature | |
| Show settings | |
| | |
| | |
| | |
| | |
| | |
| | |
| Close | Home |

Submenu "Safety controller" Select "Show safety controller settings"

and press the operating button.

The overview display shows the set-points and actual values of the main temperature controller and the safety controller and indicates the safety controller mode.

| Fixed value | | 08.03.201 | 5 05:05:06 | |
|------------------------|------|------------|------------|--|
| | | Setpoint | Actual | |
| Chamber temperature | | | 37.6 | |
| Safety controller mode | | Limit (abs | , | |
| Overtemperature | [°C] | 40.0 | 37.6 | Overview display with safety controller mode "Limit" |
| Undertemperature | [°C] | 32.0 | | (example values) |
| | | | | |
| | | | | |
| Close | | F | lome | |
| | | | | |
| Fixed value | | 08.03.201 | 5 05:05:06 | |
| | | Setpoint | Actual | |

| Fine database | | 00.00.00 | |
|------------------------|------|------------|-------------|
| Fixed value | | 08.03.20 | 15 05:05:06 |
| | | Setpoint | Actual |
| | | | |
| Chamber temperature | [°C] | 37.0 | 37.6 |
| Safety controller mode | | Offset (re | elative) |
| Overtemperature | [°C] | 3.0 | 37.6 |
| Undertemperature | [°C] | 5.0 | |
| | | | |
| | | | |
| | | | |
| Close | | | Home |

Overview display with safety controller mode "Offset" (example values)

Go back to the initial view with "Home".

18. Defrosting during refrigerating operation

BINDER cooling incubators are very diffusion-proof. To ensure high temperature precision there is no automatic cyclic defrosting device. The DCT[™] refrigerating system largely avoids icing of the evaporation plates. However, at very low temperatures the moisture in the air can condense on the evaporator plates leading to icing.



Always close the door properly.

Operation with temperature set-points above +5 °C / 41 °F at an ambient temperature of 25 °C / 77 °F:

The air defrosts the ice cover automatically. Defrosting is continually performed.

• Operation with temperature set-points below +5 °C / 41 °F:

Icing on the evaporator is possible. Defrost the chamber manually.



With temperature set-points < +5 $^{\circ}C$ / 41 $^{\circ}F$, regularly defrost the chamber manually:

• Set the temperature to 40 °C / 104 °F (in Fixed value entry mode)

Let the chamber operate for about 30 minutes with the door closed.



Excessive frosting of the evaporator is indicated by reduced refrigerating performance.

• Operation with temperature set-points below 0 °C / 32 °F

While operating the chamber with set-points below < 0 °C / 32 °F condensation is possible at the inner surface of the door around the door gasket.



In case of heavy condensation, check tightness of the door gasket.

After one or two days operation at a set-point < $0 \degree C / 32 \degree F$ a thin ice layer can cover the inner chamber door and the glass door. The amount depends of the ambient temperature and humidity. This does not influence the proper function of the refrigerating system.



Refrigerating performance decreases while operating the chamber at temperatures < 0 °C / 32 °F due to icing of the evaporators. For this reason defrost the chamber regularly, e.g. once a week.

19. Options

19.1 APT-COM[™] 4 Multi Management Software (option)

The chamber is regularly equipped with an Ethernet interface (7) that can connect the BINDER APT-COM[™] 4 Multi Management Software. The MAC Address is indicated under *Menu* > *Settings* > *Network settings* > *Show network settings* (chap. 12.9). The actual temperature, and fan speed values are given in adjustable intervals. Programming can be performed graphically via PC. Up to 100 chambers can be cross linked. For further information, refer to the APT-COM[™] 4 operating manual.

19.2 Data logger kit (option)

The BINDER Data Logger Kit offers an independent long-term measuring system for temperature, available for different temperature ranges.

The BINDER Data Logger is equipped with a keyboard and a large LCD display, alarm functions and a real-time function. Measurement data are recorded in the Data Logger and can be read out after the measurement via the RS232 interface of the Data Logger. It offers a programmable measuring interval and permits storing up to 64000 measuring values. Reading out is done with the Data Logger evaluation software. You can give out a combined alarm and status protocol directly to a serial printer.

Data Logger Kit T 220: Measuring sensor for the temperature values of the chamber: Temperature range -90 °C / 194 °F up to +220 °C / 428 °F.



For detailed information on installation and operation of the BINDER Data Logger, please refer to the mounting instructions Art. No. 7001-0204 and to the original user manual of the manufacturer, supplied with the data logger.

19.3 Zero-voltage relay alarm output (option)

With this option the chamber is equipped with a zero-voltage relay alarm contact which serves to transmit alarms to a central monitoring facility.

The connection is realized as a DIN socket. A suitable DIN plug is enclosed.

With KB / KB-UL (E4) 53, 115 the DIN (6) socket is located on the chamber rear.

With KB / KB-UL (E6) 240, 400, 720, the DIN socket (6) is located in the right lateral control panel.

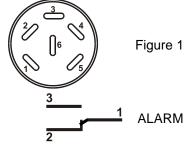



Figure 13: Pin configuration of the DIN socket (6)

Pin 1: Pole, Pin 2: Break relay, Pin 3: Make contact

In case there is no alarm, contact 1 closes with contact 3.

Closing contact 1 with contact 2 switches the zero-voltage relay alarm output.

Maximum loading capacity of the switching contacts: 24V AC/DC - 2.5A

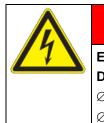

Electrical hazard through overload of contacts.

Deadly electric shock. Damage to the switching contacts and connection socket.

DANGER

 \varnothing Do NOT exceed the maximum switching load of 24V AC/DC – 2.5A.

 \varnothing Do NOT connect any devices with a higher loading capacity.

The zero-voltage relay alarm contact is activated in case of temperature tolerance range alarm and in case of a power failure.

| Displayed icon | Error when icon is displayed | Switching the alarm contact |
|----------------|-----------------------------------|-------------------------------------|
| | Power failure | immediately |
| | Temperature tolerance range alarm | 10 minutes after the error occurred |

A temperature alarm message will remain visible on the controller display during the whole time of the alarm transmission via the zero-voltage relay contact.

As soon as the cause of the alarm is rectified, you can reset the alarm transmission via the zero-voltage relay outputs together with the alarm message at the controller display by pressing the "RESET" key.

In case of a power failure, transmission of the alarm via zero-voltage relay outputs remains active for the duration of the power failure. Afterwards, both contacts will close automatically.

When using the APT-COM[™] 4 Multi Management Software (option, chap. 19.1) via the Ethernet interface of the chamber for data acquisition, the alarm message is not recorded in the APT-COM[™] protocol.

Set the tolerance limits for limit alarms by APT-COM[™] 4 separately in the AlarmCenter.

You can switch on and off the alarm output for test purpose. To access this function, select *Menu > Optional equipment > Alarm output temperature* (chap. 7.3).

19.4 Water protected internal socket (option)




The internal socket is splash proof.

IP system of protection 67 with plug or lid,100-240 V 1N \sim 50-60 Hz

Charge max. 500 W

The maximum permitted operating temperature with this option is 90 °C / 194 °F.

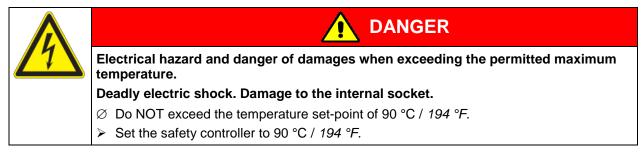




Heat emission of electrical devices connected inside the chamber may modify the temperature range.

| | NOTICE |
|----|---|
| 14 | Risk of short circuit caused by penetration of moisture into the socket. Damage to the chamber. |
| | Use the supplied plug only (IP protection type 67). Plug it in and turn it to secure contact. |
| | If the socket is not used, close the screw lid and turn it to secure. |

Turning on and off the voltage of the interior socket is possible at the chamber controller through *Menu* > *Optional equipment* > *Interior socket* >, see chap. 7.1. Turning off the incubator at the main power switch also switches the interior socket voltage-free.

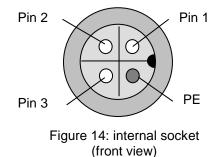

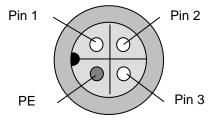


Figure 15: Supplied waterproof plug (front view)

19.5 Object temperature display with flexible Pt 100 temperature sensor (option)

You can turn on and off the object temperature display via the controller menu (chap. 7.4).

The object temperature display enables the determination of the actual temperature of the charging material during the whole process. The object temperature is measured via a flexible Pt100 temperature sensor and can be viewed at the controller display. You can immerse the sensor top protective tube of the flexible Pt 100 into liquid substances.

| Fixed value | | 08. | .03.2015 05:05:06 | |
|-------------|--------|-----------|-------------------|--------------------------------------|
| | | Setpoint | Actual value | |
| Temperature | e [°C] | 25.0 | 25.2 | |
| Fan | [%] | 100.0 | | Initial view with object temperature |
| Obj. Temp | [°C] | | 25.1 | (sample values) |
| ▶•◀ | | | | |
| User | G | uick menu | Menu | |

The object temperature data are put out together with the data of the temperature controller and can be documented by the APT-COM[™] 4 Multi Management Software (option, chap. 19.1) developed by BINDER.

Technical data of the Pt100 sensor:

- Three-wire technique
- Class B (DIN EN 60751)
- Temperature range up to 320 °C / 608 °F
- Stainless steel protective tube with a length of 45 mm / 1.78 in, material no. 1.4501

19.6 Zero-voltage relay control outputs (may be available via BINDER INDIVIDUAL Customized Solution)

The zero-voltage relay control outputs 1, 2 und 3 are used to switch any device connected via a DIN socket on the chamber rear. They permit turning on and off individually the connected devices by the controller. They can be programmed in fixed value entry mode (chap. 7.2) as well as in the time program editor (chap. 9.3.7) or the week program editor (chap. 10.3.6) via the operation lines.

The connection is realized as a DIN socket. A suitable DIN plug is enclosed.

With KB / KB-UL 53 / 115 (E4) the DIN (5) socket is located on the chamber rear.

With KB / KB-UL 240 / 400 / 720 (E6) the DIN socket (5) is located in the right lateral control panel.



Figure 16: Pin configuration of the DIN socket (5)

| Relay output 1 | Relay output 2 | Relay output 3 |
|-------------------------------|-------------------------------|-------------------------------|
| 1 Pin 1: Pin 2 Pin 2: Make | 3 Pin 3: Pin 4 Pin 4: Make | 5 Pin 5: Pin 6 Pin 6: Make |
| Switching state On: 1xx | Switching state On: x1x | Switching state On: xx1 |

Maximum loading capacity of the switching contacts: 24V AC/DC – 2.5 A



Electrical hazard through overload of contacts.

Deadly electric shock. Damage to the switching contacts and connection socket.

 \oslash Do NOT exceed the maximum switching load of 24V AC/DC – 2.5A.

 \varnothing Do NOT connect any devices with a higher loading capacity.


19.7 Interior lighting

Opening the door activates the interior lighting. When closing the door, the interior lighting goes off again.

20. Cleaning and decontamination

Clean the chamber after each use in order to prevent potential corrosion damage by ingredients of the loading material.

Prior to renewed startup, allow the chamber to completely dry after all cleaning and decontamination measures.

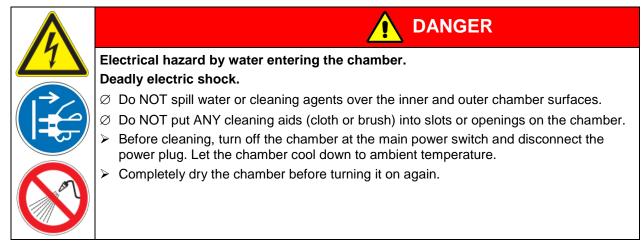

20.1 Cleaning

Disconnect the chamber from the power supply before cleaning. Disconnect the power plug.

| (A) | The interior of the chamber must be kept clean. Thoroughly remove any residues of charging material |
|-----|---|
|-----|---|

Wipe the surfaces with a moistened towel. In addition, you can use the following cleaning agents:

| Exterior surfaces inner chamber shelves door gaskets | Standard commercial cleaning detergents free from acid or halides. Alcohol based solutions. We recommend using the neutral cleaning agent Art. No. 1002-0016. | |
|---|---|--|
| Instrument panel | Standard commercial cleaning detergents free from acid or halides. We recommend using the neutral cleaning agent Art. No. 1002-0016. | |
| Zinc coated hinge parts rear chamber wall | Standard commercial cleaning detergents free from acid or halides. Do NOT use a neutral cleaning agent on zinc coated surfaces. | |

Do not use cleaning agents that may cause a hazard due to reaction with components of the device or the charging material. If there is doubt regarding the suitability of cleaning products, please contact BINDER service.

We recommend using the neutral cleaning agent Art. No. 1002-0016 for a thorough cleaning. Any corrosive damage that may arise following use of other cleaning agents is excluded from liability by BINDER GmbH. Any corrosive damage caused by a lack of cleaning, is excluded from liability by BINDER GmbH.



NOTICE

Danger of corrosion by using unsuitable cleaners. Damage to the chamber.

- $\ensuremath{\varnothing}$ Do NOT use acidic or chlorine cleaning detergents.
- \varnothing Do NOT use a neutral cleaning agent on other kind of surfaces e.g., the zinc coated hinge parts or the rear chamber wall.



For surface protection, perform cleaning as quickly as possible. After cleaning completely remove any cleaning agents from the surfaces with a moistened towel. Let the chamber dry.

Soapsuds may contain chlorides and must therefore NOT be used for cleaning.



With every cleaning method, always use adequate personal safety controls.

Following cleaning, leave the chamber door open or remove the access port plugs.



The neutral cleaning agent may cause health problems in contact with skin and if ingested. Follow the operating instructions and safety hints labeled on the bottle of the neutral cleaning agent.

Recommended precautions: To protect the eyes use sealed protective goggles. Wear gloves. Suitable protective gloves with full contact: butyl or nitrile rubber, penetration time >480 minutes.

| Danger of chemical burns through contact with skin or ingestion of the neutral cleaning agent. | |
|--|--|
| Skin and eye damage. Environmental damage. | |
| \varnothing Do not ingest the neutral cleaning agent. Keep it away from food and beverages. | |
| arnothing Do NOT empty the neutral cleaning agent into drains. | |
| Wear protective gloves and goggles. | |
| Avoid skin contact with the neutral cleaning agent. | |

20.2 Decontamination / chemical disinfection

The operator must ensure that proper decontamination is performed in case a contamination of the chamber by hazardous substances has occurred.


Disconnect the chamber from the power supply prior to chemical decontamination. Disconnect the power plug.

Do not use decontamination agents that may cause a hazard due to reaction with components of the device or the charging material. If there is doubt regarding the suitability of cleaning products, please contact BINDER service.

You can use the following disinfectants:

| Inner chamber | Standard commercial surface disinfectants free from acid or halides. |
|---------------|--|
| | Alcohol based solutions. |
| | We recommend using the disinfectant spray Art. No. 1002-0022. |

| (Jag | For chemical disinfection, we recommend using the disinfectant spray Art. No. 1002-0022. |
|------|---|
| 19 | Any corrosive damage that may arise following use of other disinfectants is excluded from liability by BINDER GmbH. |

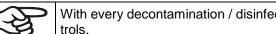

With every decontamination / disinfection method, always use adequate personal safety controls.

In case of contamination of the interior by biologically or chemically hazardous material, there are two possible procedures depending on the type of contamination and charging material.

(1) Spray the inner chamber with an appropriate disinfectant.

Before start-up, the chamber must be absolutely dry and ventilated, as explosive gases may form during the decontamination process.

(2) If necessary, have strongly contaminated inner chamber parts removed by an engineer for cleaning, or have them exchanged. Sterilize the inner chamber parts in a sterilizer or autoclave.



In case of eye contact, the disinfectant spray may cause eye damage due to chemical burns. Follow the operating instructions and safety hints labeled on the bottle of the disinfectant spray.

Recommended precautions: To protect the eyes use sealed protective goggles.





Danger of chemical burns through eye contact with the disinfectant spray.

Eye damage. Environmental damage

- \varnothing Do NOT empty the disinfectant into drains.
- Wear protective goggles.



After using the disinfectant spray, allow the chamber to dry thoroughly, and aerate it sufficiently.

21. Maintenance and service, troubleshooting, repair, testing

21.1 General information, personnel qualification

• Maintenance

See chap. 21.2.

• Simple troubleshooting

Chap. 21.3 describes troubleshooting by operating personnel. It does not require technical intervention into the chamber, nor disassembly of chamber parts.

For personnel requirements please refer to chap. 1.1.

Detailed troubleshooting

If errors cannot be identified with simple troubleshooting, further troubleshooting must be performed by BINDER Service or by BINDER qualified service partners or technicians, in accordance with the description in the Service Manual.

For personnel requirements please refer to the Service Manual.

Repair

Repair of the chamber can be performed by BINDER Service or by BINDER qualified service partners or technicians, in accordance with the description in the Service Manual.


After maintenance, the chamber must be tested prior to resuming operation.

• Electrical testing

To prevent the risk of electrical shock from the electrical equipment of the chamber, an annual repeat inspection as well as a test prior to initial startup and prior to resuming operation after maintenance or repair, are required. This test must meet the requirements of the competent public authorities. We recommend testing under DIN VDE 0701-0702:2008 in accordance with the details in the Service Manual.

For personnel requirements please refer to the Service Manual.

21.2 Maintenance intervals, service

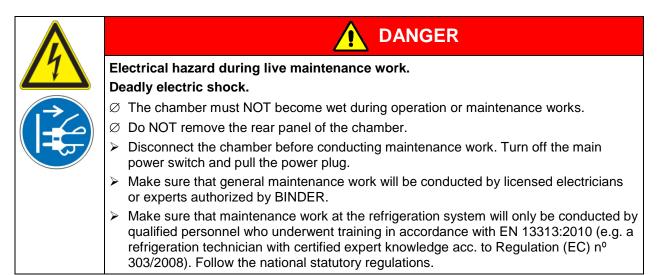

Ensure regular maintenance work is performed at least once a year and that the legal requirements are met regarding the qualifications of service personnel, scope of testing and documentation. All work on the refrigeration system (repairs, inspections) must be documented.



The warranty becomes void if maintenance work is conducted by non-authorized personnel.

BINDER

Replace the door gasket only when cold. Otherwise, the door gasket may become damaged.

With an increased amount of dust in the ambient air, clean the condenser fan (by suction or blowing) several times a year.

We recommend taking out a maintenance agreement. Please consult BINDER Service.

| BINDER telephone hotline: | +49 (0) 7462 2005 555 |
|---------------------------------------|--|
| BINDER fax hotline: | +49 (0) 7462 2005 93555 |
| BINDER e-mail hotline: | customerservice@binder-world.com |
| BINDER service hotline USA: | +1 866 885 9794 or +1 631 224 4340 x3 (toll-free in the USA) |
| BINDER service hotline Asia Pacific: | +852 390 705 04 or +852 390 705 03 |
| BINDER service hotline Russia and CIS | +7 495 988 15 16 |
| BINDER Internet website | http://www.binder-world.com |
| BINDER address | BINDER GmbH, post office box 102, D-78502 Tuttlingen |
| | |

International customers, please contact your local BINDER distributor.

21.3 Simple troubleshooting

Defects and shortcomings can compromise the operational safety of the chamber and can lead to risks and damage to equipment and persons. If there are is a technical fault or shortcoming, take the chamber out of operation and inform BINDER Service. If you are not sure whether there is a technical fault, proceed according to the following list. If you cannot clearly identify an error or there is a technical fault, please contact BINDER Service.



Only qualified service personnel authorized by BINDER must perform repair. Repaired chambers must comply with the BINDER quality standards.

| Fault description | Possible cause | Required measures |
|---------------------------|---|---|
| General | | |
| | No power supply. | Check connection to power supply. |
| Chamber without function. | | Check whether the chamber is turned on at the main power switch |
| | Wrong voltage. | Check power supply for correct voltage (chap. 4.3). |
| | Chamber fuse has responded. | Check chamber fuse and replace it if appropriate. If it responds again, contact BINDER service. |
| | Controller defective. | |
| | Nominal temperature exceeded by 10° due to chamber failure. Over temperature protective de- vice (class 1) responds. | Contact BINDER service. |



| Fault description | Possible cause | Required measures | |
|---|--|---|--|
| Heating | | | |
| Overtemperature. Chamber heating permanently, exceeding the set-point | Controller defective. | Confirm the alarm (chap. 14.4). Contact BINDER service. | |
| Notification "Heating active". | Semiconductor relay defective. | Contact BINDER Service. | |
| Alarm message "Temp. range" | Controller not adjusted. | Confirm the alarm (chap. 14.4). Calibrate and adjust controller. | |
| Excess temperature. Tolerance range alarm: Having reached the set-point, the temper- ature deviates by more than +/- 2 °C from the set-point for more than 10 minutes. Alarm message "Temp. range" | Site of installation too warm. Ambient temperature > 25 °C (chap. 3.4). | Confirm the alarm (chap. 14.4). Select cooler installation site. | |
| Excess temperature. Over- temperature safety controller class 3.1 responds. | Safety controller has turned off the heating. Limit temperature reached. Safety controller set too low. Safety controller set- point value exceeded. | Confirm the alarm (chap. 14.4). Check setting of temperature set- point and of the safety controller set-point. If appropriate, select suitable value. | |
| Alarm message: | Too much external heat load. | Confirm the alarm (chap. 14.4). Reduce heat load. | |
| "Safety controller over- | Controller defective. | | |
| temperature" | Safety controller defective. | Confirm the alarm (chap. 14.4). Contact BINDER service. | |
| | Semiconductor relay defective. | Contact BINDER Service. | |
| Excess temperature. With option safety controller class 3.3: over-temperature safety controller class 3.1 responds. | Safety controller set-point value exceeded. | Confirm the alarm (chap. 14.4). Check temperature setpoint and setting of the overtemperature safety controller. If appropriate, se- lect suitable limit value. | |
| Alarm message: | Too much external heat load. | Confirm the alarm (chap. 14.4). Reduce heat load. | |
| "Safety controller over- | Controller defective. | | |
| temperature" | Safety controller defective. | Confirm the alarm (chap. 14.4). Contact BINDER service. | |
| | Semiconductor relay defective. | | |
| Too low temperature. Chamber doesn't heat up. | Heating defective. | | |
| Notification "Heating active". | Semiconductor relay defective. | Contact BINDER service. | |
| Too low temperature. With option safety controller class 3.3: undertemperature safety controller class 3.2 responds | Safety controller set-point value exceeded. | Check setting of temperature set- point and of the safety controller class 3.2 set-point. If appropriate, select suitable value. | |
| Alarm message: | Controller defective. | Confirm the alarm (chap. 14.4). | |
| "Safety controller Under- temperature" | Safety controller defective. | Contact BINDER service. | |



| Fault description | Possible cause | Required measures | |
|--|--|--|--|
| Heating (continued) | | - | |
| range alarm: Having reached the | Chamber door not properly closed. | Confirm the alarm (chap. 14.4). Completely close chamber door. | |
| set-point, the temperature devi- ates by more than +/- 2 °C from the set-point for more than 10 minutes, or the temperature doesn't reach the tolerance range within 3 hours from turning on the chamber or closing the door. | Door gasket defective. | Confirm the alarm (chap. 14.4). Replace door gasket. | |
| Alarm message "Temp. range" | Controller defective. | Confirm the alarm (chap. 14.4). Check the controller function. If ap- propriate, contact BINDER service | |
| Too low temperature. Set-point | Chamber door not properly closed. | Completely close chamber door. | |
| temperature is not reached after specified time. | Door gasket defective. | Replace door gasket. | |
| | Controller not adjusted. | Calibrate and adjust controller. | |
| Very long heating-up times. | Chamber fully loaded. | Load the chamber less or consider longer heating-up times. | |
| Refrigerating performance | | | |
| Temperature set-point tempera- | Chamber door not properly closed. | Completely close chamber door. | |
| ture is not reached after specified time. | Door gasket defective. | Replace door gasket. | |
| | Controller not adjusted. | Calibrate and adjust controller. | |
| Low or no refrigerating | Ambient temperature above 25 °C / 77 °F (chap. 3.4). | Select cooler place of installation. | |
| performance. | Compressor not turned on. | | |
| Notification: | Solenoid valves defective. | Contact BINDER service. | |
| "Refrigeration active" | No or insufficient refrigerant. | | |
| | Too much external heat load. | Reduce heat load. | |
| Controller | 1 | | |
| Program duration longer than pro- grammed. | Inappropriate tolerances have been programmed. | For rapid transition phases, do NOT program tolerance limits in order to allow maximum heating speed. | |
| Message "Sensor rupture" | Sensor rupture between sensor and controller or Pt100 sensor defective | Contact BINDER service. | |
| No controller operation possible. Notification "Key lock activated". | Key lock activated. | Enter the key lock password (chap. 11). | |

21.4 Sending the chamber back to BINDER GmbH

If you return a BINDER product to us for repair or any other reason, we will only accept the product upon presentation of an authorization number that has previously been issued to you. An authorization number will be issued after receiving your complaint either in writing or by telephone **prior** to your sending the BINDER product back to us. The authorization number will be issued following receipt of the information below:

- BINDER product type and serial number
- Date of purchase
- Name and address of the dealer from which you bought the BINDER product
- Exact description of the defect or fault
- Complete address, contact person and availability of that person
- Exact location of the BINDER product in your facility
- A contamination clearance certificate (chap. 26) must be faxed in advance

The authorization number must be applied to the packaging in such a way that it can be easily recognized or be recorded clearly in the delivery documents.

For security reasons we cannot accept a chamber delivery if it does not carry an authorization number.

Return address:

BINDER GmbH Abteilung Service Gänsäcker 16 78502 Tuttlingen Germany

22. Disposal

22.1 Disposal of the transport packing

| Packing element | Material | Disposal |
|---|---|-------------------|
| Straps to fix packing on pallet | Plastic | Plastic recycling |
| Wooden transport box (option) | Non-wood (compressed matchwood, IPPC standard) | Wood recycling |
| with metal screws | Metal | Metal recycling |
| Pallet for KB / KB-UL 53 / 115 (E4) | Solid wood (IPPC standard) | Wood recycling |
| Pallet for KB / KB-UL 240 / 400 / 720 | Solid wood (IPPC standard) | Wood recycling |
| (E6) with foamed plastic stuffing | PE foam | Plastic recycling |
| Shipping box | Cardboard | Paper recycling |
| with metal clamps | Metal | Metal recycling |
| Top cover | Cardboard | Paper recycling |
| Removal aid for KB / KB-UL 115 | Cardboard | Paper recycling |
| | Plastic | Plastic recycling |
| Edge protection | Styropor [®] or PE foam | Plastic recycling |
| Protection of doors and racks | PE foam | Plastic recycling |
| Bag for operating manual | PE foil | Plastic recycling |
| Insulating air cushion foil (packing of optional accessories) | PE foil | Plastic recycling |

If recycling is not possible, all packing parts can also be disposed of with normal waste.

22.2 Decommissioning

Turn off the main power switch (1). Disconnect the chamber from the power supply.

When switching off the main power switch ON / OFF (1), the stored parameters remain saved.

- Temporal decommissioning: See indications for appropriate storage, chap. 3.3.
- Final decommissioning: Dispose of the chamber as described in chap. 22.3 to 22.5.

22.3 Disposal of the chamber in the Federal Republic of Germany

According to Annex I of Directive 2012/19/EU of the European Parliament and of the Council on waste electrical and electronic equipment (WEEE), BINDER devices are classified as "monitoring and control instruments" (category 9) only intended for professional use". They must not be disposed of at public collecting points.

The chambers bear the symbol for the marking of electrical and electronic equipment manufactured / placed on the market in the EC after 13 August 2005 and be disposed of in separate collection according to Directive 2012/19/EU on waste electrical and electronic equipment (WEEE) and German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG). WEEE marking: crossed-out wheeled bin with solid bar under. A significant part of the materials must be recycled in order to protect the environment.



At the end of the device's service life, have the chamber disposed of according to the German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG from 20 October 2015, BGBI. I p. 1739) or contact BINDER service who will organize taking back and disposal of the chamber according to the German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG from 20 October 2015, BGBI. I p. 1739).

NOTICE Danger of violation against existing law if not disposed of properly. Ø Do NOT dispose of BINDER devices at public collecting points. Have the device disposed of professionally at a recycling company which is certified according to the German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG from 20 October 2015, BGBI. I p. 1739).

Instruct BINDER Service to dispose of the device. The general terms of payment and delivery of BINDER GmbH apply, which were valid at the time of purchasing the chamber.

Certified companies disassemble waste (used) BINDER equipment in primary substances for recycling according to Directive 2012/19/EU. The devices must be free from toxic, infectious or radioactive substances in order to eliminate any health hazards to the employees of the recycling companies.

| (Jag | Prior to handing the chamber over to a recycling company, it is the user's responsibility that is free from toxic, infectious or radioactive substances. Prior to disposal, clean all introduced or residual toxic substances from the unit. | |
|------|---|--|
| | Prior to disposal, disinfect the unit from all sources of infection. Be aware that sources of infection may also be located outside the inner chamber. | |
| | If you cannot safely remove all toxic substances and sources of infection from the unit, dispose of it as special waste according to national law. | |
| | • Fill out the contamination clearance certificate (chap. 26) and enclose it with the unit. | |



| Danger of intoxication and infection through contamination of the chamber with toxic, infectious or radioactive substances. |
|--|
| Damages to health. |
| NEVER take a chamber contaminated with toxic substances or sources of infection for recycling according to Directive 2012/19/EU. |
| Prior to disposal, remove all toxic substances and sources of infection from the chamber. |
| A chamber from which all toxic substances or sources of infection cannot be safely re- moved must be considered as "special" waste according to national law. Dispose of it accordingly. |

The refrigerant used R134A (1,1,1,2-tetrafluorethane) is not inflammable at ambient pressure. It must not be exposed to the environment. In Europe, recovery of the refrigerant R134A (GWP 1300) is mandatory according to regulation No. 842/2006/EC. Ensure the compliance with the applicable legal requirements regarding qualification of staff, disposal, and documentation.


22.4 Disposal of the chamber in the member states of the EU except for the Federal Republic of Germany

According to Annex I of Directive 2012/19/EU of the European Parliament and of the Council on waste electrical and electronic equipment (WEEE), BINDER devices are classified as "monitoring and control instruments" (category 9) only intended for professional use". They must not be disposed of at public collecting points.

The chambers bear the symbol for the marking of electrical and electronic equipment manufactured / placed on the market in the EC after 13 August 2005 and be disposed of in separate collection according to the Directive 2012/19/EU on waste electrical and electronic equipment (WEEE). WEEE marking: crossed-out wheeled bin with solid bar under.




At the end of the device's service life, notify the distributor who sold you the device, who will take back and dispose of the chamber according to the Directive 2012/19/EU on waste electrical and electronic equipment (WEEE).

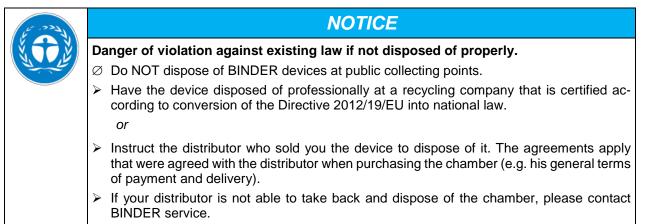

Certified companies disassemble waste (used) BINDER equipment in primary substances for recycling according to Directive 2012/19/EU. The devices must be free from toxic, infectious or radioactive substances in order to eliminate any health hazards to the employees of the recycling companies.



| (A) | Prior to handing the chamber over to a recycling company, it is the user's responsibility that it is free from toxic, infectious or radioactive substances. |
|-----|---|
| | Prior to disposal, clean all introduced or residual toxic substances from the chamber. |
| | Prior to disposal, disinfect the chamber from all sources of infection. Be aware that sources of infection may also be located outside the inner chamber. |
| | If you cannot safely remove all sources of infection and toxic substances from the chamber, dispose of it as special waste according to national law. |
| | Fill out the contamination clearance certificate (chap. 26) and enclose it with the chamber. |

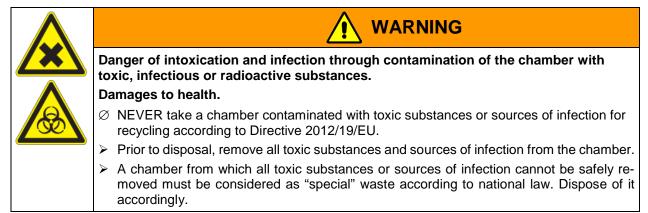

The refrigerant used R134A (1,1,1,2-tetrafluorethane) is not inflammable at ambient pressure. It must not be exposed to the environment. In Europe, recovery of the refrigerant R134A (GWP 1300) is mandatory according to regulation No. 842/2006/EC. Ensure the compliance with the applicable legal requirements regarding qualification of staff, disposal, and documentation.

22.5 Disposal of the chamber in non-member states of the EU



NOTICE

Danger of violation against existing law if not disposed of properly. Alteration of the environment.

> For final decommissioning and disposal of the chamber, please contact BINDER service.

> Follow the statutory regulations for appropriate, environmentally friendly disposal.

The main board of the cooling incubator includes a lithium cell. Please dispose of it according to national regulations.

The refrigerant used R134A (1,1,1,2-tetrafluorethane) is not inflammable at ambient pressure. It must not be exposed to the environment. In Europe, recovery of the refrigerant R134A (GWP 1300) is mandatory according to regulation No. 842/2006/EC. Ensure the compliance with the applicable legal requirements regarding qualification of staff, disposal, and documentation.

23. Technical description

23.1 Factory calibration and adjustment

This chamber was calibrated and adjusted in the factory. Calibration and adjustment were performed using standardized test instructions, according to the QM DIN EN ISO 9001 system applied by BINDER (certified since December 1996 by TÜV CERT). All test equipment used is subject to the administration of measurement and test equipment that is also constituent part of the BINDER QM DIN EN ISO 9001 systems. They are controlled and calibrated to a DKD-Standard at regular intervals.




23.2 Over current protection

The chambers are equipped with an internal fuse (over-current release) not accessible from outside. If this fuse is blown, please contact an electronic engineer or BINDER Service.


In addition, the chambers KB 53-UL / 115-UL are protected by a chamber-protection against over current, accessible from the outside. It is located at the rear of the chamber below the strain relief of the power cord. The fuse holder is equipped with a fuse clip 6.3 x32 mm. Replace this fuse only with a substitute of the same ratings. Refer to the technical data of the respective device type. If this fuse is blown, please inform an electronic engineer or BINDER service.

23.3 Definition of usable volume

The usable volume illustrated below is calculated as follows:

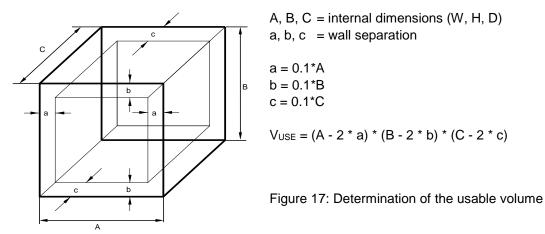

The technical data refers to the defined usable volume.

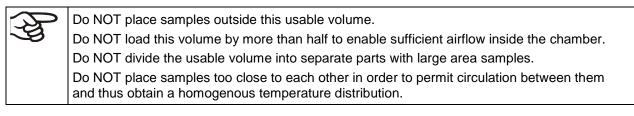

23.4 Technical data

| Chamber size | 9 | | 53 | 115 | 240 | 400 | 720 |
|---|----------------------|------------|---------------------------------|---------------------------------|--------------------------|--------------------------|---------------------------------|
| Exterior dime | | | •• | | | | |
| Width, net | | mm / inch | 635 / 25 00 | 835 / 32.88 | 925 / 36 42 | 925 / 36 42 | 1250/49.21 |
| | (incl. feet/castors) | | | 1025/40.35 | | | |
| Depth, net | | | | 650 / 25.60 | | | |
| Depth, gross | (incl. door handle, | | | 730 / 28.74 | | | |
| (/ | | mm / inch | 100 / 3.94 | 100 / 3.94 | 100 / 3.94 | 100 / 3.94 | 100 / 3.94 |
| spacers Wall clearance | e side(minimum) | mm / inch | 160 / 6.30 | 160 / 6.30 | 100 / 3.94 | 100 / 3.94 | 100 / 3.94 |
| Doors | | | 1007 0.30 | 1007 0.30 | 1007 3.34 | 1007 3.34 | 1007 3.34 |
| Number of do | ore | | 1 | 1 | 1 | 1 | 2 |
| | ner glass doors | | 1 | 1 | 1 | 1 | 2 |
| Interior dime | | [| I | I | I | I | 2 |
| Width | 11510115 | mm / inch | 100 / 15 75 | 600 / 23.62 | 650 / 25 50 | 650 / 25 50 | 070 / 28 10 |
| | | | | 480 / 18.90 | | | |
| Height | | | | 400 / 15.75 | | | |
| Depth Interior volum | • | | | | | | |
| Interior volum | | / cu.ft. | 53 / 1.9 | 115 / 4.1 | 247 / 8.7 | 400 / 14.1 | 698 / 24.6 |
| Steam space | volume | l / cu.ft. | 36 / 1.27 | 77 / 2.72 | 348 / 12.3 | 564 / 19.3 | 918 / 32.4 |
| Racks | | [| <u> </u> | | <u> </u> | <u>^</u> | |
| Quantity of racks (regular) | | | 2 | 2 | 2 | 2 | 2 |
| Quantity of racks (max.) | | 14 1 11 | 4 | 5 | 9 | 15 | 15 |
| Maximum load per rack | | Kg / Ibs | 15 / 33 | 20 / 44 | 30 / 66 | 30 / 66 | 45 / 99 |
| · · · · · · | mitted total load | Kg / Ibs | 40 / 88 | 50 / 110 | 100 / 220 | 120 / 265 | 150 / 331 |
| Weight | | F | | [| | | |
| Weight (empty | | Kg / Ibs | 78 / 172 | 106 / 234 | 171 / 377 | 221 / 487 | 304 / 670 |
| Temperature | data | | | T | | | |
| Temperature i | range | °C °F | -10 to +100 <i>14 to 212</i> | -10 to +100 <i>14 to 212</i> | -10 to +100 14 to 212 | -10 to +100 14 to 212 | -10 to +100 <i>14 to 212</i> |
| | max. | +/- K | 0,2 | 0.2 | 0,1 | 0,1 | 0,1 |
| Temperature | at 4 °C / 39.2 °F | +/- K | 0,1 | 0.1 | 0,1 | 0,1 | 0,1 |
| fluctuation | at 25 °C / 77 °F | +/- K | 0,1 | 0.1 | 0,1 | 0,1 | 0,1 |
| | at 37 °C / 98.6 °F | +/- K | 0,1 | 0.1 | 0,1 | 0,1 | 0,1 |
| | max. | +/- K | 0,1 to 1,0 | 0,1 to 1,2 | 0,1 to 0,8 | 0,1 to 1,0 | 0,1 to 0,5 |
| Temperature | at 4 °C / 39.2 °F | +/- K | 0,6 | 0.5 | 0,4 | 0,4 | 0,4 |
| uniformity (variation) | at 25 °C / 77 °F | +/- K | 0,2 | 0.1 | 0,2 | 0,2 | 0,2 |
| (variation) | at 37 °C / 98.6 °F | +/- K | 0,2 | 0.2 | 0,2 | 0,2 | 0,2 |
| Max. heat compensation at 40 °C / 104°F | | W | 150 | 210 | 300 | 500 | 500 |
| Recovery | at 4 °C / 39.2 °F | minutes | 3 | 4 | 12 | 14 | 19 |
| time after | at 25 °C / 77 °F | minutes | 1 | 1 | 1 | 1 | 1 |
| doors were open for 30 sec. | at 37 °C / 98.6 °F | minutes | 2 | 2 | 2 | 4 | 3 |

| Chamber size | | | 53 | 115 | 240 | 400 | 720 |
|--|---|-----------------|-------------------------|------------------------|-------------------------|------------------------|-------------------------|
| Electrical da | Electrical data (model versions KB053-230V, KB115-230V, KB240-230V, KB400-230V, KB720-230V) | | | | | | 20-230V) |
| IP system of | protection acc. to E | N 60529 | 20 | 20 | 20 | 20 | 20 |
| Nominal | at 50 Hz power frequency | V | 230 | 230 | 200-230 | 200-230 | 200-230 |
| voltage (+/-10%) | at 60 Hz power frequency | V | | | 200-230 | 200-230 | 200-230 |
| Current type | | | 1N~ | 1N~ | 1N~ | 1N~ | 1N~ |
| Power plug | | | shock proof plug | | | | |
| Nominal pow | /er | kW | 0.60 | 0,70 | 1.20 | 1.40 | 2.10 |
| Overvoltage | category acc. to IE | C 61010-1 | | II | II | II | II |
| Pollution deg | pree acc. to IEC 610 | 010-1 | 2 | 2 | 2 | 2 | 2 |
| Over-current B, 2 poles | release category | Amp | 16 internal | 16 internal | 16 internal | 16 internal | 16 internal |
| | ectrical data for KE ons KB053UL-120V | | | | |)V, KB720U | L-240V) |
| Nominal voltage | at 50 Hz power frequency | V | | | 100-120 | 100-120 | 200-240 |
| (+/-10%) | at 60 Hz power frequency | V | 115 | 115 | 100-120 | 100-120 | 200-240 |
| Current type | | | 1N~ | 1N~ | 1N~ | 1N~ | 2~ |
| Power plug | | NEMA | 5-15P | 5-15P | 5-20P | 5-20P | 6-20P |
| Chamber fus 250V / super | e 6,3 x 32 mm -time-lag TT | Amp | 12.5 external | 12.5 external | | | |
| Over-current B, 2 poles | Over-current release category 3, 2 poles | | 16 internal | 16 internal | 16 internal | 16 internal | 16 internal |
| Environmer KB720-230V | it-specific data (me | odel versions | KB053-230 | V, KB115-23 | 30V, KB240- | 230V, KB40 | 0-230V, |
| Noise level (| mean value.) | dB (A) | 49 | 49 | 53 | 53 | 53 |
| _ | at 4 °C / 39.2 °F | Wh/h | ≤ 260 | ≤ 250 | ≤ 350 | ≤ 380 | ≤ 420 |
| Energy con- sumption | at 25 °C / 77°F | Wh/h | ≤ 220 | ≤ 210 | ≤ 300 | ≤ 300 | ≤ 390 |
| | at 37 °C / 98.6°F | Wh/h | ≤ 70 | ≤ 75 | ≤ 270 | ≤ 330 | ≤ 360 |
| Filling weight of refrigerant R 134A (GWP 1300) | | Kg / <i>lbs</i> | 0.150 / <i>0.3</i> 3 | 0.130 / <i>0.29</i> | 0.240 / <i>0.5</i> 3 | 0.300 / <i>0.66</i> | 0.440 / <i>0.</i> 97 |
| | vironment-specificons KB053UL-120V | | | | | | L-240V) |
| Filling weigh R 134A (GW | t of refrigerant P 1300) | Kg / Ibs | 0.140 / <i>0.31</i> | 0.130 / <i>0.29</i> | 0.240 / <i>0.53</i> | 0.300 / <i>0.66</i> | 0.440 / <i>0.</i> 97 |

All technical data is specified for unloaded chambers with standard equipment at an ambient temperature of +22 °C +/- 3°C / 71.6 °F +/- 5.4 °F and a power supply voltage fluctuation of +/- 10%. Technical data is determined in accordance to BINDER Factory Standard Part 2:2015 and DIN 12880:2007.

All indications are average values, typical for chambers produced in series. We reserve the right to change technical specifications at any time.

Refrigerating performance decreases while operating the chamber at temperatures below 0 °C / 32 °F due to icing of the evaporators. For this reason defrost the chamber regularly, e.g. once a week.

If the chamber is fully loaded, the specified heating up and cooling down times may vary according to the load.

23.5 Equipment and Options (extract)

To operate the cooling incubator, use only original BINDER accessories or accessories / components from third-party suppliers authorized by BINDER. The user is responsible for any risk arising from using unauthorized accessories.

| | KB | (E4) | KB (E6) | |) |
|---|----|------|----------|----------|----------|
| Chamber size | 53 | 115 | 240 | 400 | , 720 |
| Regular equipment | | | <u> </u> | <u> </u> | |
| Multifunction program controller T4.12 with digital display | ٠ | • | • | • | • |
| Overtemperature safety controller class 3.1 acc. to DIN 12880:2007 | • | • | • | • | • |
| Ethernet interface | • | • | ٠ | • | • |
| Programmable key lock | • | ٠ | • | • | ٠ |
| Inner glass door | • | ٠ | • | • | ٠ |
| DCT™ refrigerating system with refrigerant R134a | • | ٠ | • | • | ٠ |
| Four castors (2 lockable) | | | • | • | ٠ |
| Access port 30 mm left, with silicone plug | | | • | • | • |
| Door heating | | | • | • | • |
| Options / accessories | | | | | |
| Additional shelf, stainless steel | 0 | 0 | 0 | 0 | 0 |
| Perforated shelf ,stainless steel | | | 0 | 0 | 0 |
| Reinforced rack stainless steel with 1 set of rack lockings (max. load 45 kg / <i>99.2 lb</i>) | | О | О | О | 0 |
| Reinforced rack (max. load 67 kg / 147.7 lb) with additional fixation for op- erating shakers, stirrers, roller systems | 0 | O | О | О | 0 |
| Rack lockings (4 pc.) | 0 | О | О | О | О |
| Additional access ports 10 mm / <i>0.39 in</i> / 30 mm / <i>1.18 in</i> / 50 mm / <i>1.97 in</i> or 100 mm / <i>3.94 in</i> mm with silicone plug | 0 | 0 | | | |
| Additional access ports 30 mm / <i>1.18 in</i> , 50 mm / <i>1.97 in</i> or 100 mm / <i>3.94 in</i> , with silicone plug | | | О | О | О |
| Rubber pads for safe stacking (4 pieces) | Ο | 0 | | | |
| Zero-voltage relay alarm output with DIN socket (6 poles), DIN plug in- cluded | 0 | О | О | О | О |
| Zero-voltage relay control outputs via operation lines output to DIN socket (6 poles), DIN plug included (may be available by BINDER INDIVIDUAL Customized Solutions) | 0 | o | o | О | 0 |
| Water-proof interior socket 100-240 V AC IP 67 (plug IP 67) | 0 | Ο | Ο | 0 | О |
| Object temperature display with flexible Pt 100 temperature sensor | 0 | 0 | 0 | 0 | 0 |
| Analog output for temperature 4-20mA, not adjustable, with DIN socket (6 poles), DIN plug included | 0 | 0 | 0 | 0 | 0 |
| Data Logger Kit T 220 | 0 | Ο | О | О | 0 |
| Lockable door | 0 | Ο | Ο | 0 | О |
| Over- and undertemperature safety controller class 3.3 acc. to DIN 12880:2007 | 0 | o | О | 0 | О |
| Interior lighting (KB / KB-UL 53 / 115 / 240: 15 W; KB / KB-UL 400 / 720: 30 W) | 0 | О | О | O | 0 |
| Factory calibration certificate | Ο | О | О | О | О |
| Spatial temperature measurement with certificate | 0 | 0 | О | О | О |
| Qualification folder | 0 | 0 | О | О | О |
| Stable table on wheels with castors and locking brakes | Ο | 0 | | | |

23.6 Spare parts and accessories (extract)

BINDER GmbH is responsible for the safety features of the chamber only, provided skilled electricians or qualified personnel authorized by BINDER perform all maintenance and repair, and if components relating to chamber safety are replaced in the event of failure with original spare parts. The user is responsible for any risks arising from using unauthorized accessories/components.

| | KB / KB-UL (E4) | | K | 6) | | |
|---|-----------------|-----------|-----------|-----------|-----------|--|
| Chamber size | 53 | 115 | 240 | 400 | 720 | |
| Description | ArtNo. | | | | | |
| Shelf, stainless steel | 6004-0007 | 6004-0008 | 6004-0101 | 6004-0101 | 6004-0106 | |
| Perforated shelf, stainless steel | 6004-0029 | 6004-0030 | 6004-0040 | 6004-0040 | 8009-0486 | |
| Perforated rack with add. fixation for shaker operation | 8012-0287 | 8012-0288 | 8012-0639 | 8012-0639 | 8012-0673 | |
| Reinforced rack with rack lockings | | | 8012-0638 | 8012-0638 | 8012-0674 | |
| Door gasket silicone | 6005-0095 | 6005-0096 | | | | |
| Door gasket silicone (kettle) | | | 6005-0147 | 6005-0212 | 6005-0196 | |
| Door gasket silicone (outer door) | | | 6005-0161 | 6005-0211 | 6005-0197 | |
| Intermediate door gasket silicone | | | | | 6005-0192 | |


| Description | ArtNo. |
|--|-----------|
| Rubber pads for safe stacking (4 pieces) KB / KB-UL 53 / 115 (E4) | 8012-0001 |
| Chamber fuse 6,3 x 32 mm / 250 V / 12.5 Amp super-time-lag (TT) KB 53-UL / 115-UL (E4) | |
| Securing elements for additional fastening of racks (4 pieces) KB / KB-UL 53 / 115 (E4) | 8012-0531 |
| Securing elements for additional fastening of racks (4 pieces) KB / KB-UL 240 / 400 / 720 (E6) | 8012-0620 |
| Flexible tilt protection kit | 8009-0828 |
| Plug for silicon access port d30 | 6016-0035 |
| Data Logger Kit T 220 | 8012-0715 |
| Cleaning kit (neutral cleaning agent, disinfection spray and lint-free disposable wipes, pro- tective gloves and goggles) | 8012-0503 |
| Neutral cleaning agent, 1 kg | 1002-0016 |
| Stable table on wheels with castors and locking brakes KB / KB-UL 53 / 115 (E4) | 9051-0018 |


For information on components not listed here, please contact BINDER Service.

| Validation service | ArtNo. |
|---|-----------|
| Qualification folder IQ-OQ (printed version) | 7007-0001 |
| Qualification folder IQ-OQ (digital version) | 7057-0001 |
| Qualification folder IQ-OQ-PQ (printed version) | 7007-0005 |
| Qualification folder IQ-OQ-PQ (digital version) | 7057-0005 |
| Execution of IQ-OQ | DL410200 |
| Execution of IQ-OQ-PQ | DL440500 |

| Calibration service | ArtNo. |
|---|----------|
| Calibration of temperature including certificate (1 measuring point) | DL300101 |
| Spatial temperature measurement including certificate (9 measuring points) | DL300109 |
| Spatial temperature measurement including certificate (18 measuring points) | DL300118 |
| Spatial temperature measurement including certificate (27 measuring points) | DL300127 |



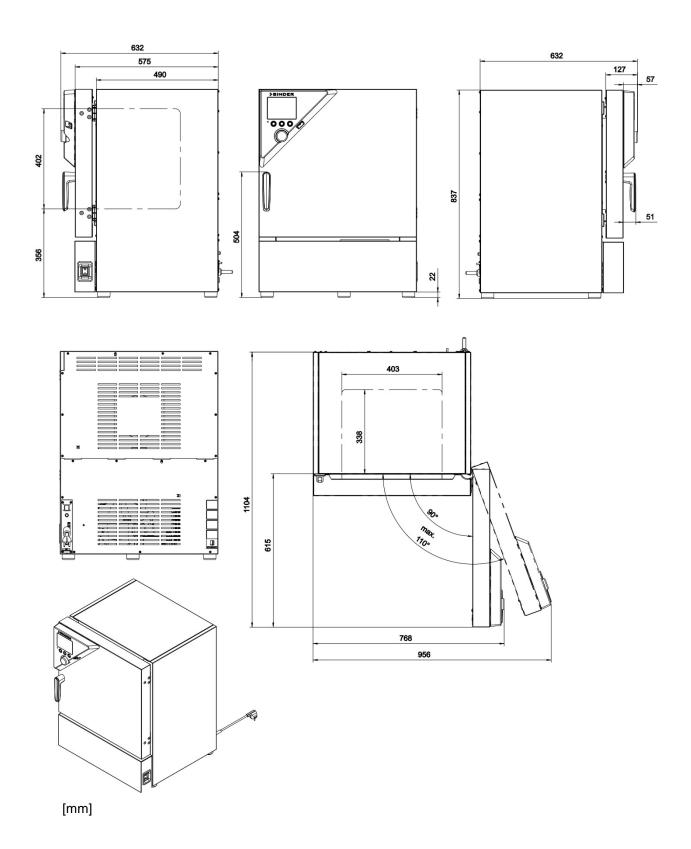


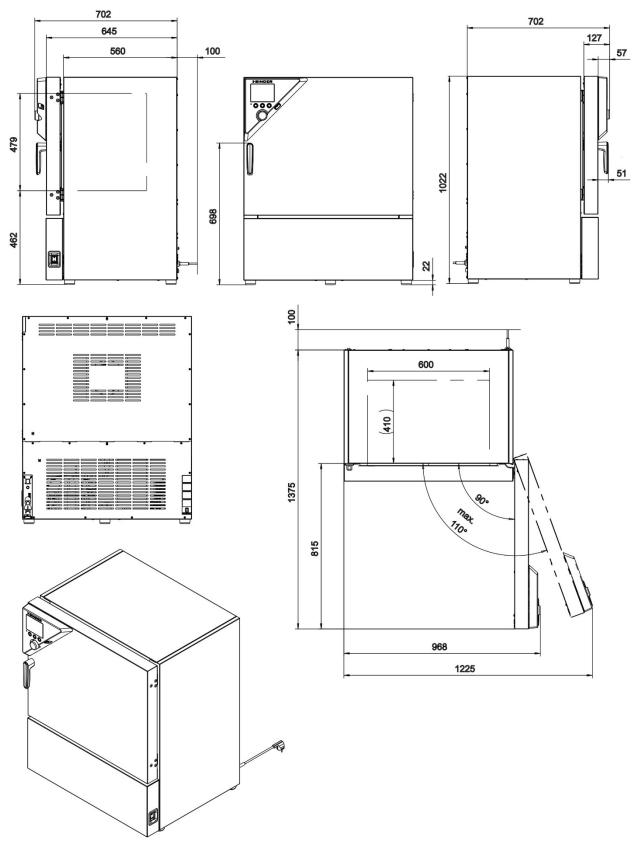




[mm]



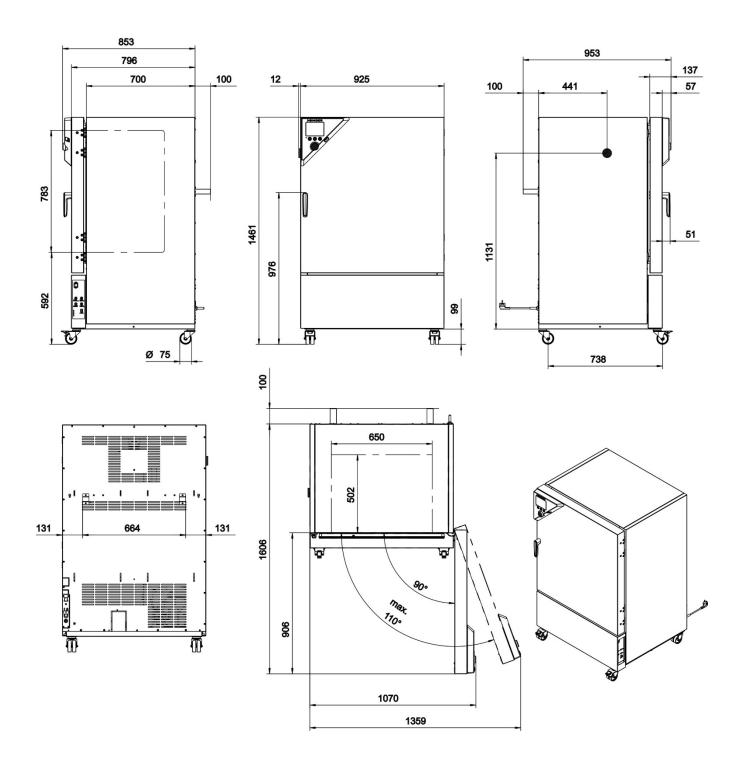


[mm]



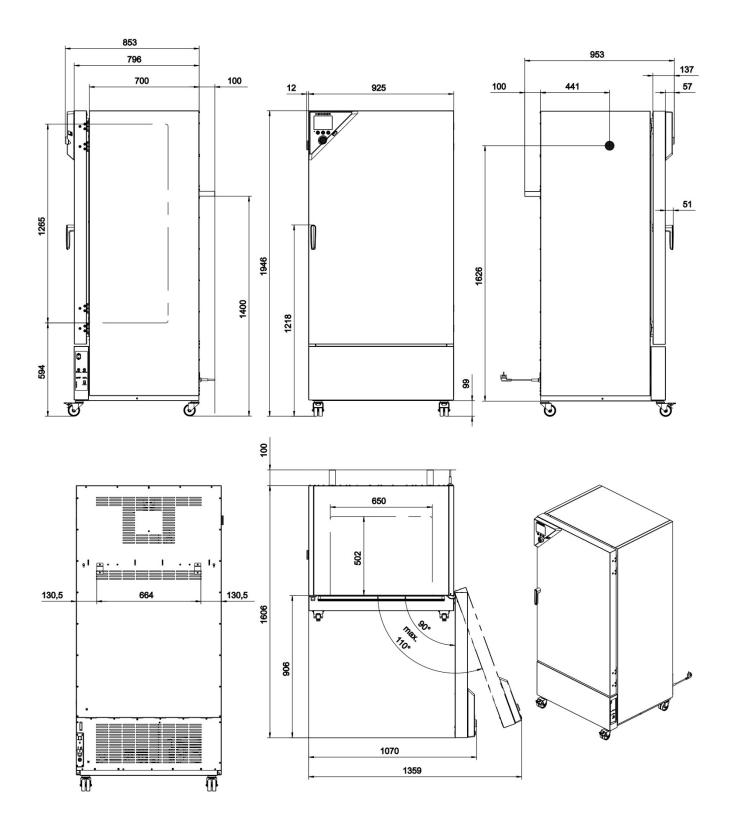

[mm]

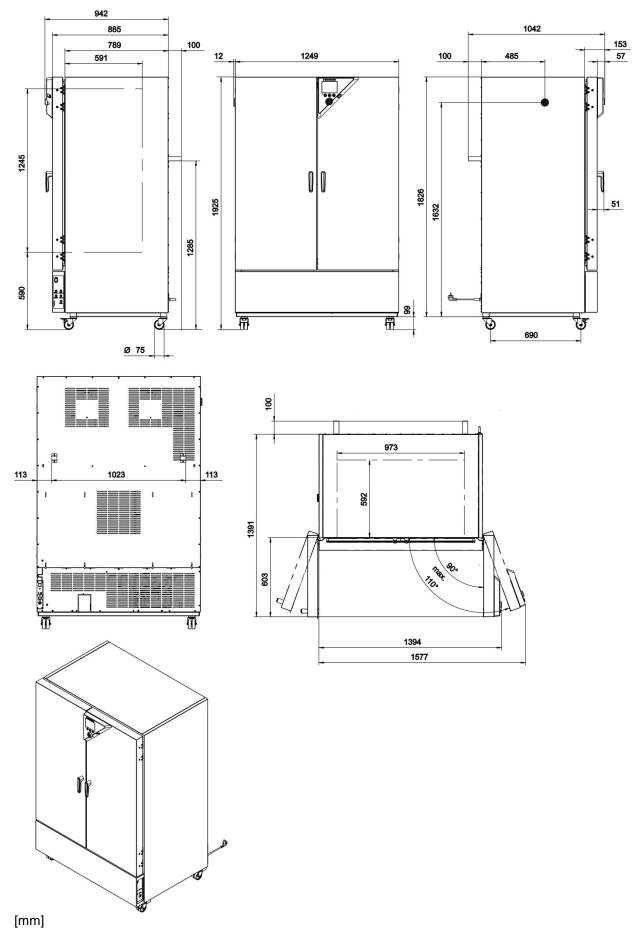

23.11 KB / KB-UL 720 dimensions

24. Certificates and declarations of conformity

24.1 EU Declaration of Conformity

| | BINDER Best conditions for your succe |
|--|--|
| СЕ EU-Konformitätserklärung / EU Ded UE / Declaración de conformidad U соответствия EU | claration of Conformity / Déclaration de conformité IE / Dichiarazione di conformità UE / Декларация |
| Hersteller / Manufacturer / Fabricant / Fabricante / Fabbricante / Производитель | BINDER GmbH |
| Anschrift / Address / Adresse / Dirección / Indirizzo / Адрес | Im Mittleren Ösch 5, 78532 Tuttlingen, Germany |
| Produkt / Product / Produit / Producto / Prodotto / Продукт | Kühlinkubatoren mit Kompressortechnologie Cooling incubators with compressor technology Incubateurs réfrigérés avec technologie de compresseur Incubadoras refrigeradas con tecnología de compressore Incubatori refrigerati con tecnologia a compressore Инкубаторы с охлаждением с компрессорной технологией |
| Typenbezeichnung / Type / Type / Tipo / Tipo / Тип | KB 53, KB 115, KB 240, KB 400, KB 720 |
| chung im Amtsblatt der europäischen Kommissior | , |
| The machines described above are in conformity Official Journal of the European Union): | with the following EC/EU Directives (as published in t |
| Les machines décrites ci-dessus sont conformes dans le Journal officiel de l'Union européenne): | s aux directives CE/UE suivantes (selon leur publicat |
| | ntes directivas de la CE/UE (publicados en el Diario ofic |
| de la Unión Europea): | |
| Gazzetta ufficiale della Commissione europea): | seguenti direttive CE/UE (secondo la pubblicazione ne |
| Машина, указанная выше, полностью с (опубликованным в Официальном журнале Евр | оответствует следующим регламентам EC/ ропейского Содружества): |
| - 2006/40/50 | ctive 2006/42/EC / Directive Machines 2006/42/EC / Direct |
| • 2014/30/EU | 2006/42/CE / Директива о машинах 2006/42/EC 014/30/EU / Directive CEM 2014/30/UE / Directiva CEM |
| Maschinenrichtlinie 2006/42/EG / Machinery directiva 2006/42/CE (Máquinas) / Direttiva macchine 2014/30/EU EMV-Richtlinie 2014/30/EU / EMC Directive 2014/30/UE / Direttiva EMC 2014/30/UE / Дире 2011/65/EU | 2006/42/CE / Директива о машинах 2006/42/EC 014/30/EU / Directive CEM 2014/30/UE / Directiva CEM ектива ЭМС 2014/30/EU Э 2011/65/EU / Directive RoHS 2011/65/UE / Directiv |
| Maschinenrichtlinie 2006/42/EG / Machinery directiva 2006/42/CE (Máquinas) / Direttiva macchine 2014/30/EU EMV-Richtlinie 2014/30/EU / EMC Directive 2014/30/UE / Direttiva EMC 2014/30/UE / Дире 2011/65/EU RoHS-Richtlinie 2011/65/EU / RoHS Directive | 2006/42/CE / Директива о машинах 2006/42/EC 014/30/EU / Directive CEM 2014/30/UE / Directiva CEM ектива ЭМС 2014/30/EU Э 2011/65/EU / Directive RoHS 2011/65/UE / Directiv Е / Директива RoHS 2011/65/EU |
| Maschinenrichtlinie 2006/42/EG / Machinery directiva 2006/42/CE (Máquinas) / Direttiva macchine 2014/30/EU EMV-Richtlinie 2014/30/EU / EMC Directive 20 (2014/30/UE / Direttiva EMC 2014/30/UE / Дире 2011/65/EU RoHS-Richtlinie 2011/65/EU / RoHS Directive | 2006/42/CE / Директива о машинах 2006/42/EC 014/30/EU / Directive CEM 2014/30/UE / Directiva CEM ектива ЭМС 2014/30/EU Э 2011/65/EU / Directive RoHS 2011/65/UE / Directiv |

| Best conditions for you | |
|---|------------------------|
| | |
| | ED |
| | |
| | |
| Die oben beschriebenen Maschinen entsprechen aufgrund ihrer Konzipierung und Bauart sowie i uns in Verkehr gebrachten Ausführung den einschlägigen grundlegenden Sicherheits- und Gesur forderungen der genannten EG/EU-Richtlinien. The machines described above are conform to the mentioned EC/EU directives in regard to the | ndheitsa |
| safety and health demands due to their conception and style of construction as well as to the ver- onto market by us. | ersion p |
| Les machines décrites ci-dessus correspondent aux demandes de sécurité et de santé des directi de la CE/UE due à leur conception et construction et dans la réalisation mise sur le marché par no | us. |
| Las máquinas descritas arriba se corresponden con los requisitos básicos pertinentes de segurida de las citadas directivas de la CE/UE debido a su concepción y fabricación, así como a la realizad da a cabo por nosotros. | ad y salı ción llev |
| Le macchine sopra descritte sono conforme ai requisiti essenziali di sanità e sicurezza pertin summenzionate direttive CE/UE in termini di progettazione, tipo di costruzione ed esecuzione mes in circolazione. | ienti de ssa da r |
| Машины описано выше, соответствует указанным директивам EC/EU в отношении тре соответствующей безопасности и здоровья по концепции и конструкции так же как и применяемая нами на рынке. | ебовані і верси |
| Die oben beschriebenen Maschinen tragen entsprechend die Kennzeichnung CE. | |
| The machines described above, corresponding to this, bear the CE-mark. | |
| Les machines décrits ci-dessus, en correspondance, portent l'indication CE. | |
| Las maquinas descritas arriba, en conformidad, llevan la indicación CE. | |
| Le macchine sopra descritte sono contrassegnate dal marchio CE. | |
| Машины описано выше, в соответствии с изложенным выше маркированы знаком СЕ. | |
| Die oben beschriebenen Maschinen sind konform mit folgenden harmonisierten Normen: | |
| The machines described above are in conformity with the following harmonized standards: | |
| Les machines décrits ci-dessus sont conformes aux normes harmonisées suivantes: | |
| Las maquinas descritas arriba cumplen con las siguientes normas: | |
| Le macchine sopra descritte sono conforme alle seguenti normative armonizzate: | |
| Машины описано выше, полностью соответствуют следующим стандартам: | |
| Sicherheit / Safety / Sécurité / Seguridad / Sicurezza / Нормативы по безопасности | |
| • EN ISO 12100:2010 + Corr. 1:2011 | |
| • EN ISO 13732-1:2008 | |
| EN 60204-1:2006 + A1:2009 + Corr. :2010 | |
| EMV / EMC / CEM / CEM / EMC / ЭМС | |
| | |
| EN 61326-1:2013 | |
| • EN 61326-1:2013 RoHS | |



BINDER Best conditions for your success

78532 Tuttlingen, 03.07.2017 BINDER GmbH

Mulinder

P. M. Binder Geschäftsführender Gesellschafter Managing Director Directeur général Director general Direttore Generale Директор

J. Bollaender

Leiter F & E und Dokumentationsbevollmächtigter Director R & D and documentation representative Chef de service R&D et autorisé de documentation Responsable I & D y representante de documentación Direttore R & D e responsabile della documentazione Глава департамента R&D представитель документации

3/3

 BINDER GmbH
 Postfach 102
 D-78502 Tuttlingen
 Address: BINDER GmbH
 Im Mittleren Ösch 5
 78532 Tuttlingen
 Germany

 Contact:
 Phone: +49 (0) 74 62 / 20 05 - 0
 |
 Fax: +49 (0) 74 62 / 20 05 - 100 |
 into@binder-world.com
 www.binder-world.com

 Managing Director:
 Dipl.-Ing. Peter M. Binder |
 District court Stuttgart, HRB 727150 |
 Company head office: Tuttlingen
 Germany

 Payment Details:
 Kreissparkasse Tuttlingen
 Account no.: 2202 611 55 |
 IBAN-Code:
 DE7464350070 0220 281155 |
 SWIFT-Code:
 SOLA DE S1TUT

 S-Account no.
 2203 611 55 |
 IBAN-Code:
 DE7464350070 0220 281155 |
 SWIFT-Code:
 DEUT DE S603

 Poutsche Bank Tuttlingen
 Account no.: 2138 709 BAN: 653 700 75 |
 IBAN-Code:
 DE56653 70075 0213870900 |
 SWIFT-Code:
 DEUT DE S603


 Recycling of old equipment according to WEEE-Reg.-no. DE 37004983
 SU
 SU
 SU
 Father School (Strateging of Strateging of Strategi



24.2 Certificate for the GS mark of conformity of the "Deutsche Gesetzliche Unfallversicherung e.V." (German Social Accident Insurance) DGUV

| Zertifikat Nr. NV 20122 vom 22.06.2020 | Prüf- und Zertifizierungsstelle Nahrungsmittel und Verpackung Fachbereich Nahrungsmittel | |
|--|--|--|
| GS-Zertifikat | | |
| Name und Anschrift des Zertifikatsinhabers: (Auftraggeber) | Binder GmbH Im Mittleren Ösch 5 78532 Tuttlingen | |
| Produktbezeichnung: | Klimaschränke Klima- und Kühlbrutschränke | |
| Тур: | KBF P 240, KBF P 720, KBF LQC 240, KBF LQC 720, KBWF 240, KBWF 720, KBF 115, KBF 240, KBF 720, KMF 115, KMF 240, KMF 720, KBW 240, KBW 400, KBW 720, KB 23, KB 53, KB 115, KB 240, KB 400, KB 720, KBF 1020, KBF 115 - DC, KBF 115 - DCL | |
| Prüfgrundlage: | GS-NV 5:2019/08 Prüfgrundsätze für Kühl- und Gefriermaschinen für Industrie und Gewerbe | |
| Zugehöriger Prüfbericht: | Prüfberichte zu den Zertifikaten NV 15127 und NV 17216 | |
| Weitere Angaben: | Das Zertifikat bezieht sich auf die im zugehörigen Prüfbericht be- schriebene Ausführung des Produkts. Nachfolgebescheinigung zu denjenigen mit den Prüfnummern NV 15127 und NV 17216 | |
| genannten Anforderungen ü GS-Zeichen an den mit dem Der Zertifikatsinhaber hat da | nmt mit den in § 21 Absatz 1 des Produktsicherheitsgesetzes berein. Der Zertifikatsinhaber ist berechtigt, das umseitig abgebildete geprüften Baumuster übereinstimmenden Produkten anzubringen. bei die umseitig aufgeführten Bedingungen zu beachten. | |
| Dieses Zertifikat einschließlie schließlich: | ch der Berechtigung zur Anbringung des GS-Zeichens ist gültig bis ein- | |
| Weiteres über die Gültigkeit, und Zertifizierungsordnung. | 21.06.2021 eine Gültigkeitsverlängerung und andere Bedingungen regelt die Prüf- | |
| | Onterschrift (Zentifizierer) | |
| Deutsche Gesetzliche Unfahlversicherung (DGUV) e / Spitzemerband der gewerblichen Berufvgemussenso | | |



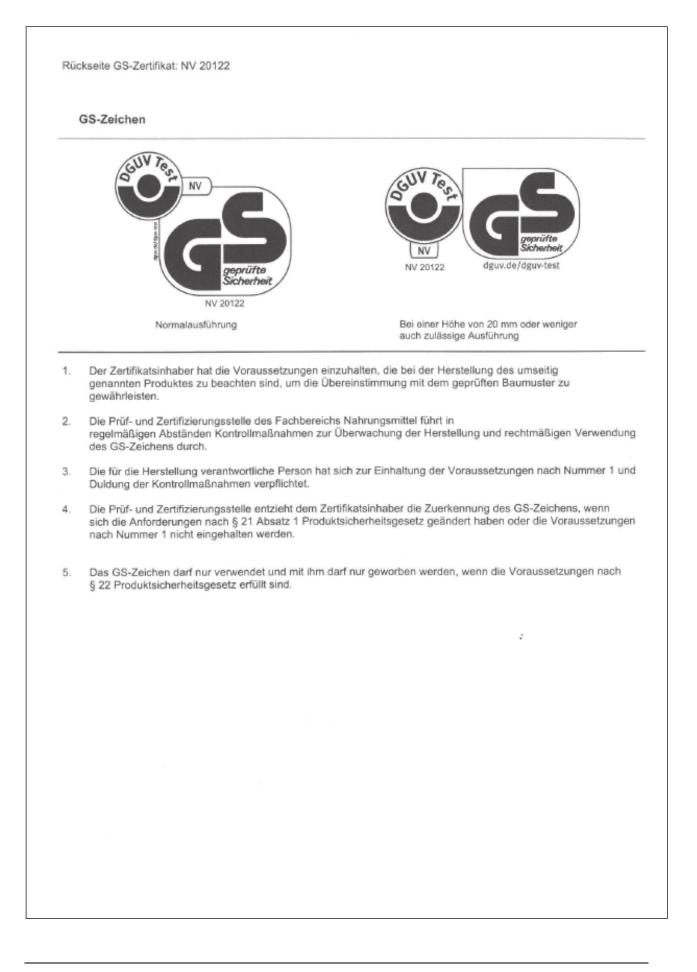

25. Product registration

Online Product Registration Register your BINDER now!

www.binder-world.com/register

The registration is free and takes just a few seconds Advantages:

- ▶ Short response times if service is needed
- ▶ Fair prices when relocating or installing equipment
- Calibration as required at no charge in case of recalls
- ► Free information on news, product upgrades and accessories

Easy registered in 3 steps:

| Nominal tampsrature XOX 1°C Not concerne XX XX Tenge, safety device XX XX Tenge, safety device XX XX You now You now You n |
|--|
| 1. List serial number here: |
| Go online: <u>www.binder-world.com/register</u> |
| 3. Register serial number |
| |
| |
| |

26. Contamination clearance certificate

26.1 For chambers located outside USA and Canada

Declaration regarding safety and health

Erklärung zur Sicherheit and gesundheitlichen Unbedenklichkeit

The German Ordinance on Hazardous Substances (GefStofV), and the regulations regarding safety at the workplace, require that this form be filled out for all products that are returned to us, so that the safety and the health of our employees can be guaranteed.

Die Sicherheit und Gesundheit unserer Mitarbeiter, die Gefahrstoffverordnung GefStofV und die Vorschriften zur Sicherheit am Arbeitsplatz machen es erforderlich, dass dieses Formblatt für alle Produkte, die an uns zurückgeschickt wird.



A completely filled out form must be transmitted via Fax (+49 (0) 7462 2005 93555) or by letter in advance, so that this information is available before the equipment/component part arrives. A second copy of this form must accompany the equipment/component part. In addition, the carrier should be informed.

Eine vollständig ausgefüllte Kopie dieses Formblattes soll per Telefax (Nr. +49 (0) 7462 2005 93555) oder Brief vorab an uns gesandt werden, so dass die Information vorliegt, bevor das Gerät/Bauteil eintrifft. Eine weitere Kopie soll dem Gerät/Bauteil beigefügt sein. Ggf. ist auch die Spedition zu informieren.

Incomplete information or non-conformity with this procedure will inevitably lead to substantial delays in
processing. Please understand the reason for this measure, which lies outside our area of influence and
will help us to speed up this procedure.

Unvollständige Angaben oder Nichteinhalten dieses Ablaufs führen zwangsläufig zu beträchtlichen Verzögerungen in der Abwicklung. Bitte haben Sie Verständnis für Maßnahmen, die außerhalb unserer Einflussmöglichkeiten liegen und helfen Sie mit, den Ablauf beschleunigen.


• Please print and fill out this form completely.

Bitte unbedingt vollständig ausfüllen!

| 1. | Unit/ component part / type: / Gerät / Bauteil / Typ: |
|-----|--|
| 2. | Serial No./ Serien-Nr.: |
| 3. | Details about utilized substances / biological substances / Einzelheiten über die eingesetzten Substanzen/biologische Materialien: |
| 3.1 | Designations / Bezeichnungen: |
| a) | |
| b) | |
| c) | |
| 3.2 | Safety measures required for handling these substances / Vorsichtsmaßnahmen beim Umgang mit diesen Stoffen: |
| a) | |
| b) | |
| c) | |
| | |

| 3.3 | Measures to be taken in case of skin contact or release into the atmosphere / Maßnahmen bei Personenkontakt oder Freisetzung: | | |
|---|--|--|--|
| a) | | | |
| b) | | | |
| c) | | | |
| d) | | | |
| 3.4 | Other important information that must be taken into account / Weitere zu beachtende und wichtige Informationen: | | |
| a) | | | |
| b) | | | |
| c) | | | |
| 4. | Declaration on the risk of these substances (please checkmark the applicable items) / Erklärung zur Gefährlichkeit der Stoffe (bitte Zutreffendes ankreuzen) : | | |
| □ 4.1 | For non toxic, non radioactive, biologically harmless materials / für nicht giftige, nicht radio- aktive, biologisch ungefährliche Stoffe: | | |
| We he Gerät/B | reby guarantee that the above-mentioned unit / component part… / Wir versichern, dass o.g. | | |
| Has not been exposed to or contains any toxic or otherwise hazardous substances / weder giftige noch sonstige gefährliche Stoffe enthält oder solche anhaften. | | | |
| | t eventually generated reaction products are non-toxic and also do not represent a hazard / auch entstandene Reaktionsprodukte weder giftig sind noch sonst eine Gefährdung darstellen. | | |
| | ntual residues of hazardous substances have been removed / evtl. Rückstände von Gefahrstoffen ent- | | |
| □ 4.2 | For toxic, radioactive, biologically harmful or hazardous substances, or any other ha- zardous materials / für giftige, radioaktive, biologisch bedenkliche bzw. gefährliche Stoffe oder an- derweitig gefährliche Stoffe. | | |
| We he | reby guarantee that … / Wir versichern, dass … | | |
| pon plet | hazardous substances, which have come into contact with the above-mentioned equipment/com- ent part, have been completely listed under item 3.1 and that all information in this regard is com- e / die gefährlichen Stoffe, die mit dem o.g. Gerät/Bauteil in Kontakt kamen, in 3.1 aufgelistet sind und alle aben vollständig sind. | | |
| | the unit /component part has not been in contact with radioactivity / das Gerät/Bauteil nicht mit Radi- ivität in Berührung kam | | |
| 5. I | Kind of transport / transporter / Transportweg/Spediteur: | | |
| Transp | ort by (means and name of transport company, etc.) Versendung durch (Name Spediteur o.ä.) | | |
| Date of | dispatch to BINDER GmbH / Tag der Absendung an BINDER GmbH: | | |
| | | | |

| We hereby declare that the following measures have been taken / Wir erklären, dass folgende Maßnah- men getroffen wurden: |
|---|
| Hazardous substances were removed from the unit including component parts, so that no hazard exists for any person in the handling or repair of these items / das Gerät/Bauteil wurde von Gefahrstoffen befreit, so dass bei Handhabung/Reparaturen für die betreffenden Person keinerlei Gefährdung besteht |
| The unit was securely packaged and properly identified / das Gerät wurde sicher verpackt und vollständig gekennzeichnet. |
| Information about the hazardousness of the shipment (if required) has been provided to the transporter / der Spediteur wurde (falls vorgeschrieben) über die Gefährlichkeit der Sendung informiert. |
| We hereby commit ourselves and guarantee that we will indemnify BINDER GmbH for all damages that are a consequence of incomplete or incorrect information provided by us, and that we will exempt BINDER GmbH from eventual damage claims by third parties./ Wir versichern, dass wir gegenüber BINDER für jeden Schaden, der durch unvollständige und unrichtige Angaben entsteht, haften und BINDER gegen eventuell entstehende Schadenansprüche Dritter freistellen. |
| We are aware that, in accordance with Article 823 of the German Civil Code (BGB), we are directly liable with regard to third parties, in this instance especially the employees of BINDER GmbH, who have been entrusted with the handling / repair of the unit / component. / Es ist uns bekannt, dass wir gegenüber Dritten – hier insbesondere mit der Handhabung/Reparatur des Geräts/des Bauteils betraute Mitarbeiter der Firma BINDER - gemäß §823 BGB direkt haften |
| |
| Name: |
| Position/Title: |
| Date / Datum: |
| Signature / Unterschrift: |
| Company stamp / Firmenstempel: |
| |
| |
| |

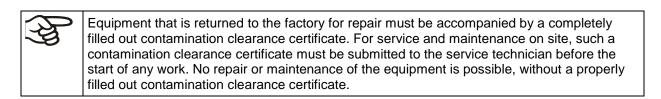

26.2 For chambers in USA and Canada

Product Return Authorization Request

Please complete this form and the Customer Decontamination Declaration (next 2 pages) and attach the required pictures. E-mail to: IDL_SalesOrderProcessing_USA@binder-world.com

After we have received and reviewed the complete information we will decide on the issue of a RMA number. Please be aware that size specifications, voltage specifications as well as performance specifications are available on the internet at <u>www.binder-world.us</u> at any time.

Take notice of shipping laws and regulations.

| | Please fill: | | |
|--------------------------------------|---------------------------|------------|---------------------------------|
| Reason for return request | O Duplicate order | | |
| | O Duplicate | e shipment | |
| | O Demo | | Page one completed by sales |
| | O Power Plug / Voltage | | 115V / 230 V / 208 V / 240V |
| | O Size does not fit space | | |
| | O Transport Damage | | Shock watch tripped? (pictures) |
| | O Other (specify below) | | |
| | | | |
| Is there a replacement PO? | O Yes | O No | |
| If yes -> PO # | | | |
| If yes -> Date PO placed | | | |
| | | | |
| Purchase order number | | | |
| BINDER model number | | | |
| BINDER serial number | | | |
| Date unit was received | | | |
| | | | |
| Was the unit unboxed? | O Yes | O No | |
| Was the unit plugged in? | O Yes | O No | |
| Was the unit in operation? | O Yes | O No | |
| Pictures of unit attached? | O Yes | O No | Pictures have to be attached! |
| Pictures of Packaging at- tached? | O Yes | O No | |

| | Customer Contact Information | Distributor Contact Information |
|---------|------------------------------|---------------------------------|
| Name | | |
| Company | | |
| Address | | |
| Phone | | |
| E-mail | | |

Customer (End User) Decontamination Declaration

Health and Hazard Safety declaration

To protect the health of our employees and the safety at the workplace, we require that this form is completed by the user for all products and parts that are returned to us. (Distributors or Service Organizations cannot sign this form)

NO RMA number will be issued without a completed form. Products or parts returned to our NY warehouse without a RMA number will be refused at the dock.

A second copy of the completed form must be attached to the outside of the shipping box.

| 1. | Unit/ component part / type: |
|-----|---|
| 2. | Serial No. |
| 3. | List any exposure to hazardous liquids, gasses or substances and radioactive material |
| 3.1 | List with MSDS sheets attached where available or needed (if there is not enough space available below, please attach a page): |
| a) | |
| b) | |
| c) | |
| 3.2 | Safety measures required for handling the list under 3.1 |
| a) | |
| b) | |
| c) | |
| 3.3 | Measures to be taken in case of skin contact or release into the atmosphere: |
| a) | |
| b) | |
| c) | |
| d) | |
| 3.4 | Other important information that must be considered: |
| a) | |
| b) | |
| c) | |

| 4. Decla | ration of Decontamination | | |
|--|--|--|--|
| For toxic, radio ous materials. | oactive, biologically and chemically harmful or hazardous substances, or any other hazard- | | |
| We hereby gu | arantee that | | |
| 4.1 Any hazardous substances, which have come into contact with the above-mentioned equipment / component part, have been completely listed under item 3.1 and that all information in this regard is complete. | | | |
| 4.2 That t | he unit /component part has not been in contact with radioactivity | | |
| | azardous substances were removed from the unit / component part, so that no hazard exists in the shipping, handling or repair of these returned unit | | |
| 4.4 The unit was securely packaged in the original undamaged packaging and properly identified on the outside of the packaging material with the unit designation, the RMA number and a copy of this declaration. | | | |
| 4.5 Shippi | ing laws and regulations have not been violated. | | |
| I hereby commit and guarantee that we will indemnify BINDER Inc. for all damages that are a con- sequence of incomplete or incorrect information provided by us, and that we will indemnify and hold harmless BINDER Inc. from eventual damage claims by third parties | | | |
| | | | |
| Name: | | | |
| Position: | | | |
| Company: | | | |
| Address: | | | |
| Phone #: | | | |
| Email: | | | |
| Date: | | | |
| Signature: | | | |

Equipment returned to the NY warehouse for repair must be accompanied by a completed customer decontamination declaration. For service and maintenance works on site, such a customer decontamination declaration must be submitted to the service technician before the start of work. No repair or maintenance of the equipment is possible without a completed form.