# ES Series Laboratory Refrigerators User's Manual

## Laboratory Freezer Range

Please read and  
keep these  
instructions

*Applicable models:*

*151F-AEV-TS*

*151F-AEW-TS*

*232F-AEV-TS*

*232F-AEW-TS*

## Contents

|  |    |
|--|----|
| Intended Use and Special Warning                       | 3  |
| Before first use                                       | 3  |
| Positioning of freezer                                 | 3  |
| Transportation and moving of freezer                   | 4  |
| Energy saving tips                                     | 4  |
| Important information                                  | 4  |
| Product information                                    | 4  |
| Product details  | 5  |
| Controller overview                                    | 6  |
| Viewing and resetting minimum and maximum temperatures | 7  |
| Changing the factory set point                         | 8  |
| Changing the temperature scale                         | 9  |
| Changing the remote alarm contacts                     | 9  |
| Alarms   | 10 |
| Remote alarm contacts                                  | 10 |
| Maintenance and Servicing                              | 12 |
| Defrosting   | 12 |
| Cleaning   | 12 |
| Reversing the door                                     | 13 |
| Storage  | 14 |
| Electrical Connection                                  | 15 |
| Disposal   | 16 |
| Troubleshooting  | 17 |
| Warranty   | 18 |
| Contact Information                                    | 18 |

## Intended Use and Special Warning

This equipment is tested in compliance with established regulations and then shipped ready for use. This equipment is intended for general purpose cold storage and is not to be used in the presence of explosive gases or mixtures and / or close to sources with high magnetic or electric fields. Not intended for flammable material storage.

*If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.*

Thank you for purchasing this ES series cold storage equipment. Before you first use this laboratory freezer please carry out the following actions:

- Check that the freezer has not been damaged in any way during transportation.
- If any damage is found it must be reported your local dealer immediately.
- Ensure all packaging has been removed, including cardboard, polystyrene and any tape used to hold shelves in place for transportation.
- The freezer has been cleaned prior to despatch; however, we advise that it should be cleaned using lukewarm water containing a mild detergent and a soft cloth.
- We recommend that the freezer should be left in an upright position for 24 hours prior to switching on.
- This equipment is not intended for use by persons (including children) with reduced physical, sensory or mental capabilities, or lack of experience and knowledge, unless they have been given supervision or instruction concerning use of the equipment by a person responsible for their safety.
- Children should not be allowed to play with the equipment.

### Positioning and safely using your ES series Freezer

- **Warning** – to reduce risk of the unit falling over, do not tip any unit beyond 10 degrees
- Make sure the equipment is placed in a dry, well ventilated site, away from heat sources.
- The equipment must be placed on a level surface. If necessary, adjust the feet on the product so the equipment remains level during operation.
- The equipment should have sufficient clearance around it to provide adequate ventilation – 10cm at each side, 6cm at the rear and 2.5cm at the top.
- Before ANY maintenance activity, pull the plug out of the socket by gripping the plug, DO NOT pull on the plug cord.
- DO NOT use double-wire extension cords. If an extension cord is necessary make sure it is a cord, which possess a protection safety certificate.
- If the main power supply cord is damaged, it should be replaced by a qualified service engineer.
- DO NOT attempt any repairs to the equipment yourself (apart from those highlighted in the Maintenance and Service Section). Repairs carried out by someone without the relevant training are putting their personal safety at risk and will invalidate your warranty.
- Do not store explosive substances such as aerosol cans with a flammable propellant in this equipment.
- **WARNING:** this equipment is only to be operated by authorised personnel or personnel who have correct knowledge on how the equipment operates.

## Transporting and moving your ES series Freezer

The freezer must always be moved in the vertical position. The cabinet must not be tilted anymore than 10°. In the event that the cabinet is tilted more than 10°, the power supply must not be connected until the equipment has stood upright for at least 24 hours.

### Energy saving tips

This equipment is cooled by energy-efficient R600a refrigerant. To keep running costs as low as possible always:

- Position the equipment away from heat sources.
- Make sure the air can circulate freely around the equipment, don't block ventilation grid.
- Ensure that products being stored are below room temperature upon entry.
- Make sure the door is opened as little as possible whilst in use and closed as quickly as possible to prevent unnecessary temperature fluctuations.

### Important Information

The freezer contains environment-friendly, non ozone depleting refrigerant R600a. As R600a is a flammable gas, it is important to avoid damage to the refrigeration circuit during transport and installation. If the refrigeration circuit is damaged, avoid using a naked flame in the vicinity of the freezer and connecting power to the freezer. Also make sure there is good ventilation in the room. If you are in doubt please contact your supplier.

- **WARNING:** Keep ventilation openings, in the equipment enclosure or in the built-in structure, clear of obstruction.
- **WARNING:** Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended by the manufacturer.
- **WARNING:** Do not damage the refrigerant circuit
- **WARNING:** Do not use electrical equipments inside the food storage compartments of the equipment, unless they are of the type recommended by the manufacturer.

### Product Information

This freezer has been designed and built with the intention of the product being used within a Laboratory and is designed to store contents between -12 and -20°C.

Please see below for specific product information:

| Model | Height (mm) | Width (mm) | Depth (mm) | Weight (Kg) | Refrigerant type | Ambient/Room working temperature | Door Type |
|-------|-------------|------------|------------|-------------|------------------|----------------------------------|-----------|
| 151F  | 845         | 595        | 635        | 36          | R600a            | +10°C to +32°C                   | Solid     |
| 232F  | 1565        | 595        | 670        | 60          |                  |                                  | Solid     |

Product details

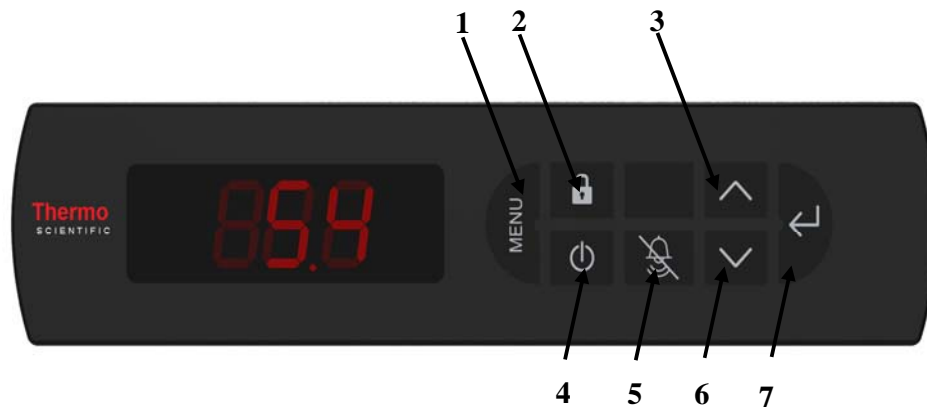
1. Lock
2. Electronic Controller
3. Door Handle
4. Wire Shelf
5. Wire Basket
6. Adjustable foot (2 at front of equipment)

There are also 2 roller castors at the rear of the equipment to ease product movement.

Please contact your local representative for more information on replacement parts

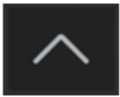
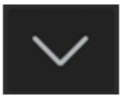
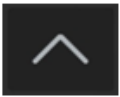
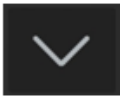
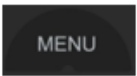
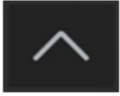
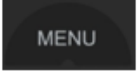
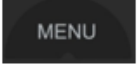
## Controller overview and operation

1. Menu
2. Lock keypad
3. Maximum temperature / Up
4. Power
5. Mute
6. Minimum temperature / Down
7. Set / Enter

## Viewing & resetting minimum and maximum temperatures

The factory default maximum temperature is XXX °C and the minimum temperature is XXX °C.

| Function   | Press   | Hold button(s) down for | Example screen  |
|--|---|-------------------------|---|
| To view Maximum air temperature<br><i>N.B. Displays maximum since last reset</i> |    | 3 seconds               |    |
| To view Minimum air temperature<br><i>N.B. Displays minimum since last reset</i> |    | 3 seconds               |    |
| To reset air temperature<br><i>N.B. This will delete all previous values</i>     |       | 3 seconds               |    |
| To view Maximum "load" temperature   |    | 3 seconds               |    |
| followed by  |    | 1 second                |    |
| followed by<br><i>N.B. Displays maximum "load" temperature since last reset</i>  |   | 3 seconds               |  |
| To view Minimum "load" temperature   |    | 3 seconds               |  |
| followed by  |    | 1 second                |  |
| followed by<br><i>N.B. Displays minimum "load" temperature since last reset</i>  |    | 3 seconds               |  |
| To reset "load" temperature<br><i>N.B. This will delete all previous values</i>  |    | 3 seconds               |  |
| followed by  |    | 1 second                |  |
| followed by  |   | 3 seconds               |  |

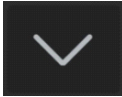
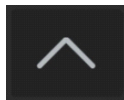
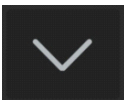
### Changing the factory set point

(the set point is the point at which the freezer normally holds the temperature at; in rare circumstances such as high/low room temperatures it may be necessary to adjust this slightly. Changing the factory setting is not recommended)

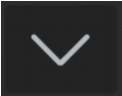
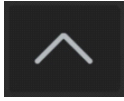
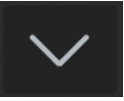
| Function   | Press  | Hold button(s) down for | Example screen   |
|--|--|-------------------------|--|
| To change the set point  |   | 3 seconds               |                         |
|  | followed by   | Once                    |                         |
|  | followed by   | 1 second                | <br>(current set point) |
|  | followed by  / <br>(Increase / Decrease) | up to 10 times          |                        |
|  | followed by   | 1 second                |                       |
| Note: set point can be increased or decreased in +/- 0.1° increments by a MAXIMUM of 1°C |  |                         |  |

### Changing the temperature scale between Celsius (°C) and Fahrenheit (°F)

| Function                    | Press   | Hold button(s) down for | Example screen  |
|-----------------------------|---|-------------------------|---|
| To change temperature scale |    | 3 seconds               |                                    |
| followed by                 |    | Two times               |                                    |
| followed by                 |    | 1 second                | <br><small>(current scale)</small> |
| followed by                 |  /  | Once                    |                                    |
| followed by                 |    | 1 second                |                                    |

### Changing the remote alarm contacts

| Function                          | Press   | Hold button(s) down for | Example screen  |
|-----------------------------------|---|-------------------------|---|
| To change the alarm contact state |    | 3 seconds               |                                    |
| followed by                       |    | Three times             |                                    |
| followed by                       |    | 1 second                | <br><small>(current state)</small> |
| followed by                       |  /  | Once                    |                                    |
| followed by                       |    | 1 second                |                                    |

Note: NO = Normally Open, NC = Normally Closed

## Warning Alarms

This freezer has been designed to provide both an audible and visual alarm should any of the following situations occur:

- Door is left open for a set period
- Internal air temperature goes outside the set parameters, either too high or too low
- Internal 'load' temperature goes outside the set parameters, either too high or too low

Should one of these alarms occur this can be muted by pressing  on the controller.

The 'load' temperature is a simulated temperature which is intended to provide an indication of the temperature of the freezer's contents (e.g. a 5ml sample). The temperature of a freezer's contents will change more slowly (up or down) than the air temperature within that freezer should there be a fault or the door is left open.

## Remote alarm contacts

All ES series laboratory freezers include remote alarm contacts and these terminal contacts enable the user to connect to a remote control station; these contacts can be found on the rear of the cabinet. **WARNING:** These contacts must only be accessed by a competent electrician.

In order to use these contacts the following must be carried out:

1. The freezer **MUST** be disconnected from the power supply before removing the terminal cover.
2. Unscrew the cover (figure 1) on the right hand side of the rear of the cabinet to expose the connectors (figure 2).

Figure 1

Figure 2

3. Use the right hand outer positions that correspond to the left hand cable positions for your connections using 0.75mm cable stripped to 0.6mm.
4. Depress the larger slot above the cable entrance on the right hand side with a small electrical screwdriver. This allows the cable to be inserted into the slot directly below.

In the event of a door open alarm, high or low temperature alarm or power failure a remote relay will switch within the controller to make an electrical circuit.

***Please note that the contacts supplied are voltage free. A voltage of between 12V DC and 230V AC can be connected at the contacts. The maximum load must not exceed 2A. The minimum power rating is 500mA / 12V AC.***

In a normal condition (i.e. no alarm) the relay OPENS the contact for the remote alarm facility.  
In an alarm condition the relay CLOSES the contact for the remote alarm facility.

## Maintenance and Service of your ES Series Freezer

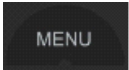
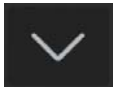
### Defrosting

All laboratory freezers require manual. In certain conditions (e.g. when the freezer is very full or when any new items are stored in the freezer) frost may form on the rear wall as the compressor must run for long periods of time.

Before manually defrosting the cabinet, make sure the power lead is disconnected from the mains. Any loose frost found on the evaporator can be removed carefully using a wooden or plastic scrapper.

**WARNING:** Do not use any knives or sharp metallic objects to remove frost.

Alternatively a non-scheduled defrost can sometimes be carried out as follows:

| Function                             | Press   | Hold button(s) down for | Example screen  |
|--------------------------------------|---|-------------------------|---|
| To carry out a non-scheduled defrost |    | 3 seconds               |    |
| followed by                          |   | Four times              |   |
| followed by                          |  | 1 second                |  |
| followed by                          |  | 1 second                |  |

N.B. This does not replace the automatic defrost function on relevant models

### Cleaning

- Your freezer should be cleaned regularly.
- Before cleaning you **MUST** switch off and disconnect your freezer from the power supply.
- The whole freezer except the door gasket can be cleaned with mild cleaning detergent.
- The door gasket should be cleaned with water only and wiped dry.
- The handle should be cleaned regularly with soap and water.
- Remove shelves/basket/door furniture and wipe with a soft cloth. **DO NOT** put them in a dishwasher.
- Detergents containing abrasives or acids are not suitable for the cleaning.
- When all cleaning has been completed and accessories replaced, reconnect the plug with **dry hands**.

## Reversing the door

In order to reverse the door please follow the below instructions. This should be carried out by 2 people for safety purposes.

1. Tilt the freezer backwards by 45° and lean on a vertical, hard surface such as a wall.
2. Remove the feet from the bottom, right hand hinge (figures 1 and 2).
3. Remove the screws from the hinge (figure 3) and remove hinge.

Figure 1

Figure 2

Figure 3

4. Pull the door from the bottom to remove from the freezer.
5. Remove the top, right hand hinge (figure 4).
6. Install the hinge on the top, left hand side (figure 5).

Figure 4

Figure 5

7. Remove the bottom, left hand feet and base plate (figures 6 and 7).

Figure 6

Figure 7

8. Insert the door into the top, left hand hinge (figure 8).
9. Install the bottom, left hand hinge (figure 9) and screw in the feet.

Figure 8

Figure 9

10. Add the feet and base plate to the bottom, right hand side.
11. Remove the caps from the screws holding the handle in position.
12. Remove the handle screws.
13. Remove the caps on the opposite side of the door.
14. Screw the handle into position and add the caps to cover the screws.

### Storage

If the freezer is to remain unused for a period of time (e.g. during refurbishment work) switch the equipment off by pressing the power button  and disconnect the power supply to the cabinet at the wall socket. It is also advised that the equipment is unplugged from the power supply. Empty, defrost and clean the freezer. Make sure the door is left slightly ajar to prevent any bad odours forming within the cabinet.

## Electrical Connection

The equipment is intended for connection to alternating current. The connection values for voltage (V) and frequency (Hz) are declared on the name plate in the cabinet. Power must be connected via a wall socket with a switch. It is strongly recommended that the wall socket is easily accessible. All earthing requirements stipulated by the local electricity authority. The cabinet plug and wall socket should give correct earthing. If in doubt, contact your local supplier or an authorised electrician.

If the supply cord is damaged, it must be replaced by the manufacturer, its service agent or similarly qualified persons in order to avoid a hazard.

### **WARNING: THIS EQUIPMENT MUST BE EARTHED**

The flexible cord (mains lead) fitted to this equipment has three cores for use with a 3-pin 13 amp plug. If a BS 1363 (13-amp) fused plug is used it should be fitted with a 13-amp fuse.

The cores in the mains lead are coloured in accordance with the following code:

**GREEN AND YELLOW: EARTH**

**BLUE: NEUTRAL**

**BROWN: LIVE**

These colours might not correspond with the colour markings identifying the terminals in your plug.

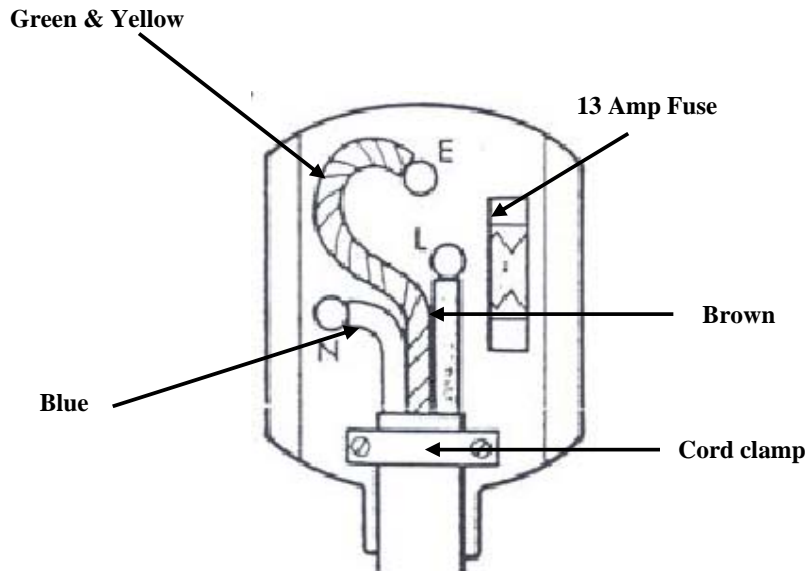
Note:

First Connection to the mains

The equipment provides information about the higher temperature (Hi alarm) till the moment of cooling.

The alarm is signalled by a sound signal and “Hi” caption on the display. To cancel the alarm, press any button.

Until the moment of cooling, the alarm is repeated every 15 minutes. Such an operating mode is normal until the required temperature settings are reached.

### Disposal of your freezer

When disposing your refrigerator do not take it to regular municipal waste containers. Instead, take it to an electrical/electronic waste recycling and re-use centre. This refrigerator contains insulation gases and refrigerant and must be disposed of in the appropriate manner. Please contact the manufacturer or local authority for advice on how best to dispose of this product.

A relevant label has been placed on the refrigerators packaging (see below). The product has been manufactured of recyclable materials. Ask your local environmental care authority for information concerning location of such facilities.

## Troubleshooting

| <b>PROBLEM</b>  | <b>REASON</b>  | <b>SOLUTION</b>  |
|---|--|--|
| The equipment does not work   | A gap in the electric system circuit   | <ol style="list-style-type: none"> <li>1. Make sure the plug is properly inserted into the socket</li> <li>2. Make sure the socket is not faulty</li> <li>3. Make sure the power supply cord is not damaged</li> </ol> |
|   | The compressor hardly ever activates   | <ol style="list-style-type: none"> <li>1. Check to see if the ambient temperature is below +10C</li> </ol>   |
| The temperature inside the equipment is not low enough (the compressor operates continuously) | The door doesn't shut tight or is opened too often   | <ol style="list-style-type: none"> <li>1. Rearrange the products so they don't hamper the door</li> <li>2. Shorten the time the door is open for</li> </ol>  |
|   | The ambient temperature is above that which has been outlined in the product specification                 | <ol style="list-style-type: none"> <li>1. Check that the equipment is operating at an ambient temperature that meets the product specification</li> </ol>  |
|   | Air circulation at the back of the equipment is hampered   | <ol style="list-style-type: none"> <li>1. Move equipment further away from the wall</li> </ol>   |
|   | The equipment is placed such that it is subject to constant sunlight, or another heat source e.g. Radiator | <ol style="list-style-type: none"> <li>1. Move the equipment to a different place away from heat sources or direct sunlight</li> </ol>   |
| Water collects at the bottom part of the refrigerator   | The contents touch the back wall of the cabinet  | <ol style="list-style-type: none"> <li>1. Move the products so they aren't in contact with the back wall</li> </ol>  |
|   | The drain opening is clogged   | <ol style="list-style-type: none"> <li>1. Take the cleaning plug and unclog the condensate opening</li> </ol>  |
| The equipment makes too much noise  | The equipment is not level   | <ol style="list-style-type: none"> <li>1. Place the equipment on an even surface or use adjustable feet</li> </ol>   |
|   | The equipment is touching another object   | <ol style="list-style-type: none"> <li>1. Reposition the equipment so it is detached from any other objects</li> </ol>   |

## WARRANTY

### **Thermo Scientific ES Series Refrigerators and Freezers European Warranty Policy**

---

---

#### **Thermo Scientific ES Series Refrigerators • European Warranty • 12 Months Full Warranty Parts and Labor**

During the first twelve (12) months from shipment, Thermo Fisher Scientific Inc, through its authorized Dealer or service organizations, will at its option and expense repair or replace any part found to be non-conforming in material or workmanship. Thermo Fisher Scientific Inc reserves the right to use replacement parts, which are used or reconditioned. Replacement or repaired parts will be warranted for only the unexpired portion of the original warranty.

This warranty does not apply to damage caused by (i) accident, misuse, fire, flood or acts of God; (ii) failure to properly install, operate or maintain the products in accordance with the printed instructions provided, (iii) causes external to the products such as, but not limited to, power failure or electrical power surges, (iv) improper storage and handling of the products, (v) use of the products in combination with equipment or software not supplied by Thermo Fisher; or (vi) installation, maintenance, repair, service, relocation or alteration of the products by any person other than Thermo Fisher or its authorized representative. To obtain proper warranty service, you must contact the nearest authorized service center or Dealer. Thermo Fisher Scientific, Inc's own shipping records showing date of shipment shall be conclusive in establishing the warranty period. At Thermo Fisher's option, all non-conforming parts must be returned to Thermo Fisher postage paid and replacement parts are shipped FOB Thermo Fisher's location.

#### **Limitation of Liability**

THIS WARRANTY IS EXCLUSIVE AND IN LIEU OF ALL OTHER WARRANTIES, WHETHER WRITTEN, ORAL, OR IMPLIED. NO WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE SHALL APPLY. THERMO FISHER DOES NOT WARRANT THAT THE PRODUCTS ARE ERROR-FREE OR WILL ACCOMPLISH ANY PARTICULAR RESULT. THERMO FISHER SHALL NOT BE LIABLE FOR ANY INDIRECT OR CONSEQUENTIAL DAMAGES INCLUDING, WITHOUT LIMITATION, DAMAGES TO LOST PROFITS OR LOSS OF PRODUCTS.

## CONTACT INFORMATION

**For technical or sales support please contact your local distributor or Thermo Scientific at the following numbers:**

**Europe:** Austria +43 1 801 40 0, Belgium +32 2 482 30 30, France +33 2 2803 2180, Germany national toll free 08001-536 376,

Germany international +49 6184 90 6940, Italy +39 02 02 95059 434-254-375, Netherlands +31 76 571 4440, Nordic/Baltic countries +358 9 329 100,

Russia/CIS +7 (812) 703 42 15, Spain/Portugal +34 93 223 09 18, Switzerland +41 44 454 12 12, UK/Ireland +44 870 609 9203

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