
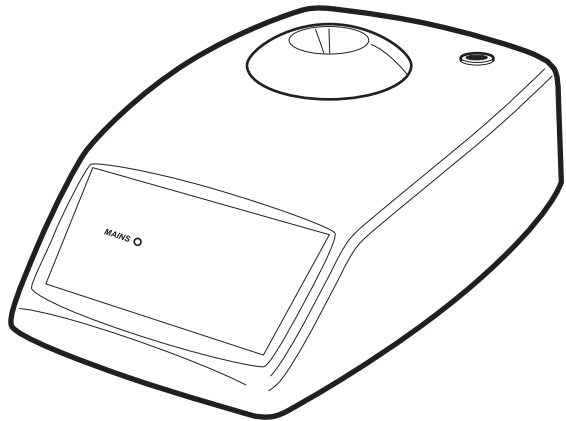
# Cole-Parmer®

## V-100 / V-200 Series

### Vortex Mixers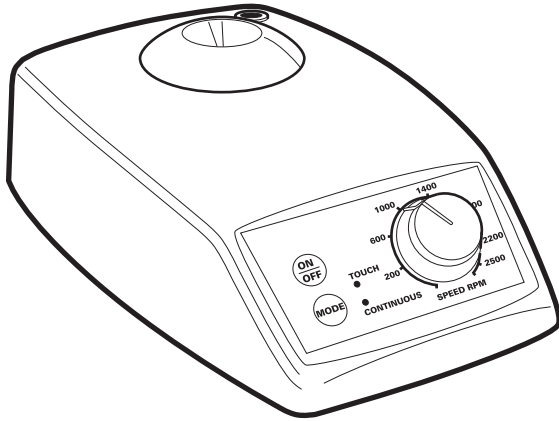

Instruction Manual  
SA7022 Version 1.4

**Cole-Parmer®**  
essentials

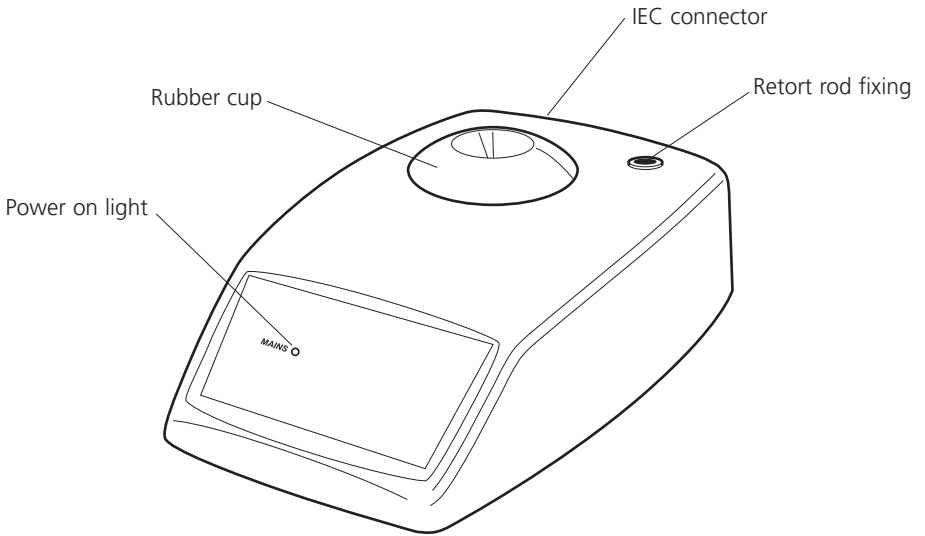
Figure 1: V-100 Fixed Speed model

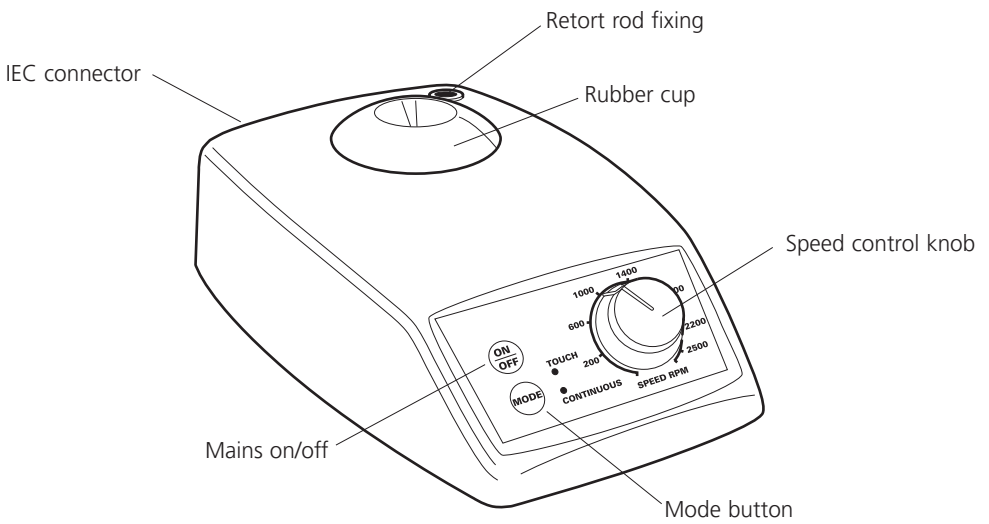
Figure 2: V-200 Variable Speed model

Thank you for purchasing this Cole-Parmer product. To get the best performance from the equipment, and for your personal safety, please read these instructions carefully before use.

This equipment is designed to operate under the following conditions:-

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature range +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuations not exceeding 10%
- ❖ Overvoltage category II IEC60364-4-443
- ❖ Pollution degree 2 IEC664
- ❖ Use with a minimum distance all round of 200mm from walls or other items

If the equipment is not used in the manner described in this manual and with accessories other than those recommended by the manufacturer, the protection provided may be impaired.

## Electrical Installation

**THIS EQUIPMENT MUST BE EARTHED**

Before connection please ensure that the line supply corresponds to that stated on the rating label.

Power requirements:

Model	Wattage
V-100	20W
V-200	20W

The Vortex Mixer is provided with a NEMA 5-15 plug for 120 V installations and a UK 3-pin and a "Schuko" 2-pin plug for 230 V installations.. Choose the lead appropriate for your electrical installation and discard the other.

Should neither lead be suitable, take the lead with the U.K. plug and replace the plug with a suitable alternative.

**THIS OPERATION SHOULD ONLY BE UNDERTAKEN BY A QUALIFIED ELECTRICIAN**

**NOTE:** Refer to the equipment's rating plate to ensure that the plug and fusing are suitable for the voltage and wattage stated.

The wires in the power cable (230 V) are coloured as follows:

BROWN	- HOT/LIVE
BLUE	- NEUTRAL
GREEN/YELLOW	- EARTH

The wires in the power cable (120 V) are coloured as follows

BLACK	- HOT/LIVE
WHITE	- NEUTRAL
GREEN	- EARTH

**IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN**

Should the mains lead need replacing, a cable of 1mm<sup>2</sup> of harmonised code H05W-F connected to an IEC320 plug should be used.

The appropriate mains lead should be connected BEFORE connection to the mains supply.

### Safety Advice Before Use

- ❖ Never move or carry the unit when in use or connected to the mains electricity supply.
- ❖ DO NOT position the product so that it is difficult to access the ON/OFF switch.
- ❖ DO NOT position the product so that it is difficult to disconnect it from the mains supply using the mains plug.
- ❖ The mains outlet socket used should be located close to the equipment and readily identifiable and accessible to users.
- ❖ Do not mix flammable liquids or use the equipment in hazardous atmospheres.
- ❖ In the case of mains interrupting, a fault or mechanical failure, the unit will NOT continue to operate on removal of fault.
- ❖ Mechanical energy can lead to breakage of glass vessels. Use with care.

## Operation

Position the apparatus on a firm level surface, ensuring that the IEC connector is in easy reach.

V-100 Fixed speed version (see Fig 1)

Connect the unit to the mains. Green neon lights. To mix tube, press base firmly down onto the rubber cap, the unit will start. Once the contents of tube are mixed, lift clear, the unit will stop until another tube is pressed onto the cup.

Description	Catalogue Number
Microtitre plate holder	SA8/1
Spare rubber cup	SA7013
Retort rod, 300mm	SR2

### V-200 Variable speed version (see fig 2)

Ensure that the speed control knob is fully to the left. Connect the unit to the mains.

Turn ON by pressing the <On / Off> button once (a green neon lights to show the unit is in the touch mode).

To mix tube, press base firmly down onto the rubber cap. Once the contents of tube are mixed, lift clear. Select the optimum mixing speed by turning the control knob to the right.

To select continuous mode, press the <mode> button (the light will change to show continuous mode). The unit will now remain on until either the <mode> button is pressed to select touch mode or the unit is switched OFF.

### Accessory heads

N.B. Accessory heads are for use with model V-200 only. Accessory heads should always be used in the continuous mode.

To change the head, remove the rubber cup by pulling upwards and store in a safe place. Insert the accessory head ensuring that the two slots in the head are in line with the lugs in the vortex mixer top.

The microtitre plate holder (SA8/1) should not be used at speeds above 800rpm.

For a comprehensive list of parts required by service engineers conducting internal repairs and a service manual, please contact the Technical Service Department of Antylia Scientific quoting both the model and serial number.

Only spare parts supplied or specified by Antylia Scientific or its agents should be used. Fitting of non-approved parts may affect the performance and safety features designed into the instrument.

If in any doubt, please contact the Technical Service Department of Antylia Scientific or the point of sale.

Antylia Scientific Ltd.  
Beacon Road,  
Stone, Staffordshire,  
ST15 0SA, United Kingdom  
Tel: +44 (0)1785 812121  
Email: cpservice@antylia.com  
Web: www.coleparmer.com

## Maintenance & Servicing

**WARNING:** Ensure the unit is disconnected from the mains electricity supply before attempting maintenance or servicing.

Periodically clean the instrument using a damp cloth and mild detergent solution. Do not use harsh or abrasive cleaning agents.

Any repairs or replacement of parts **MUST** be undertaken by suitably qualified personnel.

The following spares and accessories are available from your laboratory dealer.

# Technical Specification

	V-100	V-200
Speed, rpm	2,500	250-2,500
Orbit, mm	4.2	4.2
Continuous mode	No	Yes
Accessory heads	No	Yes
Maximum load	1 Kg	
Overall dimension	78 x 135 x 215mm	
Net weight	3.2 kg	
Electrical supply	90 - 240V, 20W	

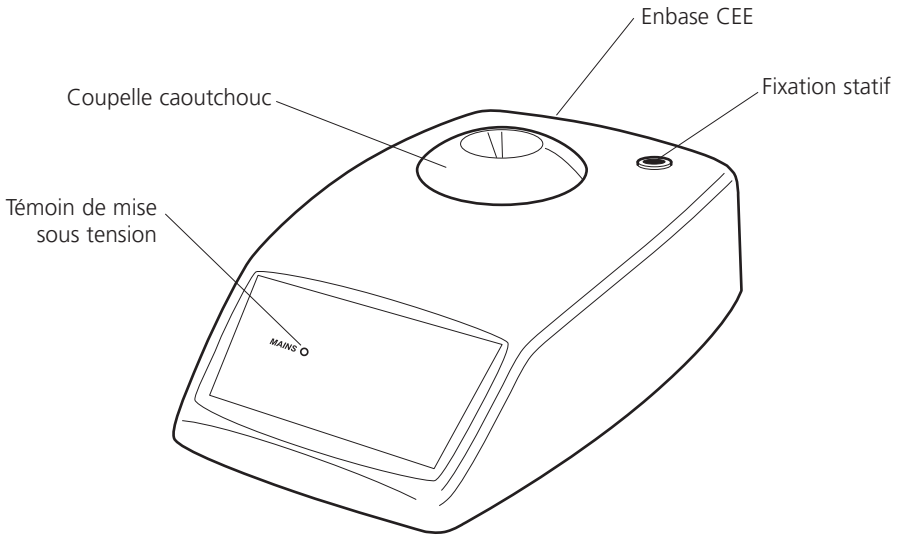
Figure 1: Version vitesse constante

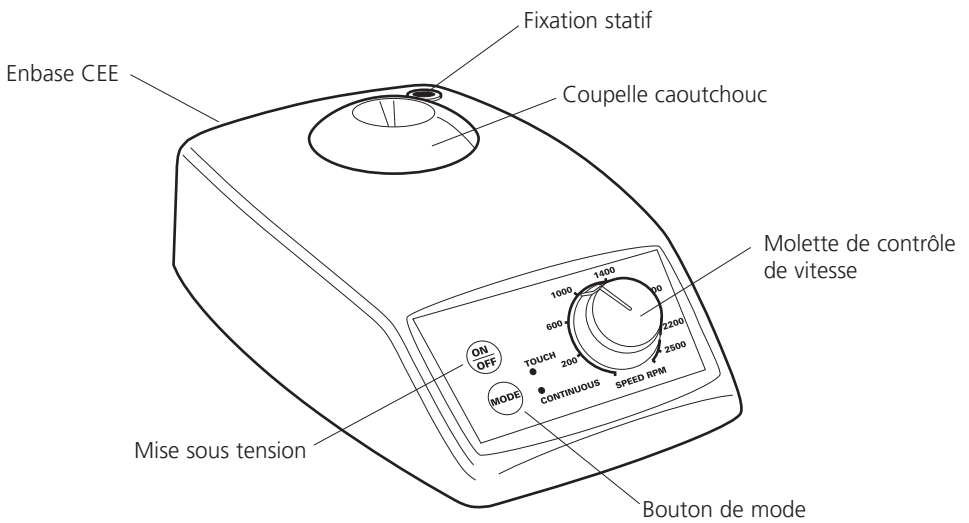
Figure 2: Version vitesse variable

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# Notes

**CE UK CA** This product meets the applicable CE Directives and UKCA Legislation for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in its vicinity will meet these standards and so we cannot guarantee

that interference will not occur in practise. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advise before use, contact the manufacturer.

Declaration of Conformity is available to view online at [www.coleparmer.com](http://www.coleparmer.com)

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T: +1 847 549 7600

**Cole-Parmer®**  
essentials

**Antylia Scientific Ltd.**  
Beacon Road,  
Stone,  
Staffordshire,  
ST15 0SA,  
United Kingdom