*Atomic Spectrometry*

# ***AAnalyst 800***

***Atomic Absorption Spectrometer***

***User's Guide***

***PERKIN ELMER***

Bodenseewerk Perkin-Elmer GmbH is registered for the design and manufacture of laboratory analytical equipment under the quality requirements of BS EN ISO 9001.

Certificate No. FM 22178

### **Notice**

The information contained in this document is subject to change without notice.

Software programs are protected by copyright. It is unlawful to duplicate these programs in any manner other than for the user's personal use. Specifically, it is unlawful to use or duplicate these programs other than for use with the purchaser's original computer.

### **Release history**

<b>Part Number</b>	<b>Release</b>	<b>Publication Date</b>
0993-5256	1	February 1998
	2	May 1998
	3	September 1998
	4	June 1999

### **Trademarks**

Perkin Elmer is a registered trademark of affiliates of Perkin Elmer LLC.

AA WinLab, AAnalyst, Lumina, and THGA are trademarks of affiliates of Perkin Elmer LLC.

Windows is a trademark and Microsoft is a registered trademark of Microsoft Corporation.

Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

### **Copyright information**

This document contains proprietary information that is protected by copyright.

All rights are reserved. No part of this document may be reproduced in any form whatsoever or translated into any language without the prior written permission of the Perkin-Elmer Corporation or one of its subsidiaries.

Copyright ©1998–1999 affiliates of Perkin Elmer LLC

Printed in the Federal Republic of Germany

Technical Documentation

Bodenseewerk Perkin-Elmer GmbH

D-88647 Ueberlingen, Federal Republic of Germany

# Contents

---

	<i>page</i>
<b>Chapter 1 Regulatory Information</b>	
Customer Service .....	1-3
Company Name and Addresses .....	1-3
Supplies, replacement parts, and accessories .....	1-3
Regulatory information .....	1-4
IEC 1010 compliance .....	1-4
Electrical protection .....	1-4
Electromagnetic compatibility (EMC) .....	1-5
Safety practices and conventions .....	1-6
Correct use of the instrument .....	1-7
Electrical safety .....	1-8
Environment .....	1-10
Symbols used on the instrument .....	1-11
Warning markings on the instrument .....	1-12
Warning markings at the front of the spectrometer .....	1-12
Warning markings on the rear panel .....	1-14
Warning markings on the deuterium lamp cover .....	1-16
<b>Chapter 2 Safety Information</b>	
Introduction .....	2-3
Correct use of analytical instruments .....	2-3
Laboratory hygiene .....	2-4
Laboratory ventilation .....	2-4
Safe handling of chemicals .....	2-5
Waste disposal .....	2-7
Contents of waste containers .....	2-7
Disposing of defective lamps .....	2-7
UV radiation .....	2-8
0993-5256	<b>C-1</b>

	<b>page</b>
Magnetic field .....	2-8
Zeeman graphite furnace systems .....	2-8
High temperatures .....	2-9
Burner system .....	2-9
Graphite furnace .....	2-9
Quartz tube atomizer cell .....	2-9
Handling compressed gases .....	2-10
Summary of gas hazards .....	2-10
Identifying cylinders .....	2-10
Storing cylinders .....	2-11
Handling cylinders .....	2-12
Safety practices for flame atomization .....	2-13
Safety interlocks .....	2-13
Safe use of burner gases .....	2-13
Air supply .....	2-14
Nitrous oxide .....	2-14
Acetylene .....	2-15
Safe operation of the flame .....	2-16
The drain system .....	2-17
Flashbacks .....	2-17
Hazards with flame atomization .....	2-18
Safety practices for electrothermal atomization .....	2-19
Safety interlocks .....	2-19
Inert gas .....	2-19
Safety checks .....	2-19
Hazards with electrothermal atomization .....	2-20
Safety practices for FIAS and mercury/hydride systems .....	2-21
FIAS-Cells – To all users who have an MHS-20 in addition to FIAS .....	2-21
Safety checks .....	2-21
Hazards with FIAS and mercury/hydride systems .....	2-22
References for laboratory safety practice .....	2-23

	<i>page</i>
<b>Chapter 3 Spectrometer Installation</b>	
Before you install the spectrometer . . . . .	3-3
Moving and reinstalling the spectrometer . . . . .	3-4
Preparing the spectrometer to be moved . . . . .	3-4
Moving the spectrometer . . . . .	3-5
Reinstalling the spectrometer . . . . .	3-5
Connecting the burner gases . . . . .	3-6
Connecting the furnace gases . . . . .	3-9
Connecting the drain system to the burner . . . . .	3-14
Installing the furnace autosampler . . . . .	3-14
Installing the cooling system . . . . .	3-15
Electrical connections . . . . .	3-18
<b>Chapter 4 Preparing the System for Analyses</b>	
Switching on the system . . . . .	4-3
Installing lamps . . . . .	4-7
Installing sample trays . . . . .	4-11
Shutting down the system . . . . .	4-12
Emergency shutdown . . . . .	4-14
<b>Chapter 5 Spectrometer: System Maintenance</b>	
Perkin Elmer Service . . . . .	5-3
Troubleshooting . . . . .	5-3
Maintenance checklist . . . . .	5-4
Spectrometer maintenance procedures . . . . .	5-5
Cleaning the windows of the optical system . . . . .	5-5
Maintaining the electrical components . . . . .	5-7
Changing the air filters . . . . .	5-9
Replacing the deuterium lamp . . . . .	5-10
<b>Chapter 6 Parts and Supplies</b>	
Parts provided . . . . .	6-3
Replacement parts . . . . .	6-5
Obtaining supplies, replacement parts, and accessories . . . . .	6-5

	<i>page</i>
<b>Chapter 7 Laboratory Requirements</b>	
Introduction .....	7-3
Space requirements for the spectrometer system .....	7-4
Magnetic field .....	7-7
Electrical requirements .....	7-8
Environmental conditions .....	7-9
Fume ventilation .....	7-10
Gas supply requirements .....	7-11
Burner gases .....	7-12
Furnace gases .....	7-14
Cooling the spectrometer system .....	7-15
Lamps .....	7-16
Computer and printer .....	7-16
<b>Chapter 8 Spectrometer: System Description</b>	
The AAnalyst 800 Spectrometer System .....	8-3
The optical system .....	8-6
The gas and cooling connectors .....	8-8
The electrical connectors and switches .....	8-10
Technical data .....	8-13
<b>Translations</b>	
Symbols used on the instrument .....	T-3
Warning messages .....	T-11
<b>Index</b>	

<b>Contents</b>	<b>page</b>
Customer Service .....	1-3
Company Name and Addresses .....	1-3
Supplies, replacement parts, and accessories .....	1-3
Regulatory information .....	1-4
IEC 1010 compliance .....	1-4
Electrical protection .....	1-4
Electromagnetic compatibility (EMC) .....	1-5
Safety practices and conventions .....	1-6
Correct use of the instrument .....	1-7
Electrical safety .....	1-8
Environment .....	1-10
Symbols used on the instrument .....	1-11
Warning markings on the instrument .....	1-12
Warning markings at the front of the spectrometer .....	1-12
Warning markings on the rear panel .....	1-14
Warning markings on the deuterium lamp cover .....	1-16

## Customer Service

### Company Name and Addresses

**U.S.A.:**

Perkin Elmer LLC  
761 Main Avenue  
Norwalk, Connecticut 06859-0010  
U.S.A.  
Phone: 800-762-4000 or  
(+1) 203-762-4000  
Fax: (+1) 203-762-4228

**Germany:**

EG&G Perkin Elmer  
Verkauf und Service  
Rengoldshauser Strasse 11  
D-88662 Ueberlingen  
Germany  
Phone: (49-7551) 919-0  
Fax: (49-7551) 919-149

**United Kingdom:**

Perkin Elmer Ltd.  
Post Office Lane  
Beaconsfield, Bucks, HP9 1QA  
England  
Phone: (44-1494) 676161  
Fax: (44-1494) 67-9331 & -9333

**Perkin Elmer on the Internet:**

<http://www.perkin-elmer.com/ai>

### Supplies, replacement parts, and accessories

Supplies, replacement parts, and accessories can be ordered directly from Perkin Elmer, using the part numbers quoted in the guides provided with the instrument.

Perkin Elmer's catalog service offers a full selection of high-quality supplies.

To place an order for supplies and many replacement parts, request a free catalog, or ask for information:

- If you are located *within the U.S.*, call toll free **1-800-762-4002**, 8 a.m. to 8 p.m. EST. Your order will be shipped promptly, usually within 24 hours.
- If you are located *outside of the U.S.*, call your local Perkin Elmer sales or service office.

## **Regulatory information**

This instrument has been designed to comply with a wide variety of international standards governing the safety of laboratory equipment. In routine use, the instrument poses virtually no risk to you. If you take some simple, common-sense precautions, you can maintain the continued safe operation of the instrument.

### **IEC 1010 compliance**

This instrument has been designed and tested in accordance with EN 61010-1:1993 (IEC 1010-1): *Safety requirements for electrical equipment for measurement, control, and laboratory use*, including Amendments 1 and 2 to this standard, and EN 61010-2-061:1996 (IEC 1010-2-061:1995): *Particular requirements for laboratory atomic spectrometers with thermal atomization and ionization*.

### **Electrical protection**

#### **Insulation**

Class I as defined in IEC 1010-1.

#### **Installation Category**

This instrument is able to withstand transient overvoltage according to Installation Category II as defined in IEC 1010-1 and IEC 664.

#### **Pollution Degree**

This equipment will operate safely in environments that contain nonconductive foreign matter and condensation up to Pollution Degree 2 as defined in IEC 1010-1 and IEC 664.

## **Electromagnetic compatibility (EMC)**

### **European Union (EMC Directives)**

This instrument has been designed and tested to meet the requirements of the EU Directives 89/336/EEC and 92/31/EEC. It complies with the generic EMC standard EN 61326-1:1997 *Electrical equipment for measurement, control, and laboratory use – EMC requirements – General requirements*.

This instrument has passed the following EMC tests:

<i>Emission of conducted and radiated noise</i>	<i>Electromagnetic compatibility</i>
EN 55011:96	EN61000-4-2:95
EN 61000-3-2:95	EN61000-4-3:95
EN 61000-3-3:95	EN61000-4-4:95
	EN61000-4-5:95
	EN61000-4-6:96

### **United States (FCC)**

This instrument is classified as a digital device used exclusively as industrial, commercial, or medical equipment. It is exempt from the technical standards specified in Part 15 of FCC Rules and Regulations, based on Section 15.103 (c).

### **Japan (FCC)**

This instrument has been tested and found to comply with the limits of a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense. Changes or modifications not expressly approved by the manufacturer could void the user's authority to operate the equipment.

## Safety practices and conventions

The guides provided with the instrument contain information and warnings that must be followed by the user to ensure safe operation and to maintain the instrument in a safe condition. This advice is intended to supplement, not supersede, the normal safety code of behavior prevailing in the country of operation.

The information provided does not cover every safety procedure that should be practiced. Ultimately, maintenance of a safe laboratory environment is the responsibility of the user and the user's organization.

Further safety practices are described in the chapter 'Safety Information'.

Possible hazards that could harm the user or result in damage to the instrument are clearly stated at appropriate places throughout these guides.

Any of the following safety conventions can be used throughout these guides:

This symbol alerts you to situations that could result in **personal injury** to yourself or other persons.  
Details about these circumstances are in a box like this one.

This symbol alerts you to the risk of **electric shock** that could result in **personal injury** to yourself or other persons.  
Details about these circumstances are in a box like this one.

This symbol alerts you to the risk of **hot surfaces** that could cause **personal injury** to yourself or other persons.  
Details about these circumstances are in a box like this one.

This symbol alerts you to the risk of **ultraviolet radiation** that could cause **eye damage** to yourself or other persons.  
Details about these circumstances are in a box like this one.

---

**Caution:**

The term **CAUTION** alerts you to situations that could result in **serious damage to the instrument** or other equipment.

Details about these circumstances are described in a message similar to this one.

---

### **Correct use of the instrument**

Before you install or use your instrument, and in order to get the best results, you should be familiar with all of the instruments in the system and know how to operate them. You should also be aware of the safety procedures in force in your laboratory, especially those concerning atomic spectroscopy instruments. Read the guides supplied with the instruments before you start.

If you use the instrument in a manner not specified in the guides, or if you use it for a purpose other than that intended, you may damage the instrument, or compromise your own, or someone else's, safety.

This instrument should only be operated by persons who are suitably qualified and have received adequate training.

## **Electrical safety**

To ensure satisfactory and safe operation of the instrument, it is essential that the green/yellow lead of the line power cord is connected to true electrical earth (ground).

### **Warning: Electrical Hazard**

W2.1

*Any interruption of the protective conductor inside or outside the instrument or disconnection of the protective conductor (earth/ground) terminal is likely to make the instrument dangerous.*

*Intentional interruption is prohibited.*

### **Lethal voltages are present in the instrument**

- *Even with the power switch off, line power voltages can still be present within the instrument.*
- *When the instrument is connected to line power, terminals may be live, and opening covers or removing parts (except those to which access can be gained without use of a tool) is likely to expose live parts.*
- *Capacitors inside the instrument may still be charged even if the instrument has been disconnected from all voltage sources.*

**Warning: Unauthorized Adjustments and Servicing**

W1.2

Only a Perkin Elmer service engineer or similarly trained and authorized person should be permitted to service the instrument.

- Do not attempt to make adjustments, replacements, repairs, or modifications to this instrument except as described in the documentation supplied with the instrument.

When working with the instrument:

- Connect the instrument to a correctly installed line power outlet that has a protective conductor (earth/ground).
- Do not attempt to make internal adjustments or replacements except as directed in the guides provided with the instrument.
- Do not operate the instrument with any covers or parts removed.
- Disconnect the instrument from all voltage sources before opening it for any adjustment, replacement, maintenance, or repair.  
If, afterwards, the opened instrument must be operated for further adjustment, maintenance, or repair, this must only be done by a skilled person who is aware of the hazard involved.
- Use only fuses with the required current rating and of the specified type for replacement.  
**Do not use makeshift fuses or short-circuit the fuse holders.**
- Whenever it is likely that the instrument is no longer electrically safe for use, make the instrument inoperative and secure it against any unauthorized or unintentional operation.

The instrument is likely to be electrically unsafe when it:

- Shows visible damage.
- Fails to perform the intended measurement.
- Has been subjected to prolonged storage under unfavorable conditions.
- Has been subjected to severe transport stresses.

## **Environment**

### **Operating conditions**

**Warning: Explosive Atmosphere**

W1.3

*This instrument is not designed for operation in an explosive atmosphere.*

***The instrument will operate correctly under the following conditions:***

- Indoors.
- Ambient temperature +15 °C to +35 °C (+59 °F to +95 °F).
- Ambient relative humidity 20% to 80%, without condensation.
- Altitude in the range 0 m to 2000 m.

***You can store the instrument safely under the following conditions:***

- Ambient temperature –20 °C to +60 °C (–4 °F to +140 °F).
- Ambient relative humidity 20% to 80%, without condensation.
- Altitude in the range 0 m to 2000 m.

When you remove the instrument from storage and before you put it into operation, allow it to stand for at least a day under the approved operating conditions.

## Symbols used on the instrument

Any of the following symbols can be used on the instrument:

This symbol indicates that there is a potential hazard and that the user must look in the guide for an explanation of the hazard and how to avoid it.

This symbol indicates the risk of electric shock.

This symbol indicates the risk of hot surfaces.

This symbol indicates the off position of the main power switch or circuit breaker. With the switch in this position, the instrument is disconnected entirely from the line power supply.

This symbol indicates the on position of the main power switch or circuit breaker. With the switch in this position, the instrument is connected to the line power supply but is not necessarily switched on and operational.

This symbol indicates the operational on/off switch. This switch turns the instrument on and off, but it does not disconnect the instrument from the line power supply.

This symbol indicates alternating current.

This symbol indicates that there is a potential hazard due to a strong magnetic field and that persons wearing heart pacemakers should remain at least 0.6 m away from the furnace.

## Warning markings on the instrument

### Warning markings at the front of the spectrometer

**A**

*Risk of hot surfaces*

*Risque de surfaces chaudes*

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*Oberflächen*

**B**

**Warning: Moving Parts – Risk of Injury**

*Some moving parts of the instrument are accessible in normal operation.*

*Keep hands, clothing and other objects away from the moving parts of the instrument.*

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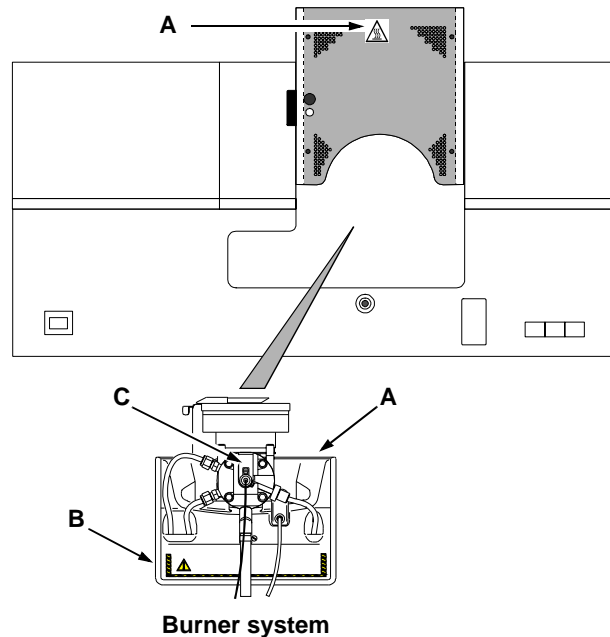
**C**

*Nebulizer clamp must be lowered and locked over the nebulizer flange.*

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**B**

**Warning: Moving Parts – Risk of Injury**

*Some moving parts of the instrument are accessible in normal operation.*  
*Keep hands, clothing and other objects away from the moving parts of the instrument.*

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**D**

**Strong magnetic field**

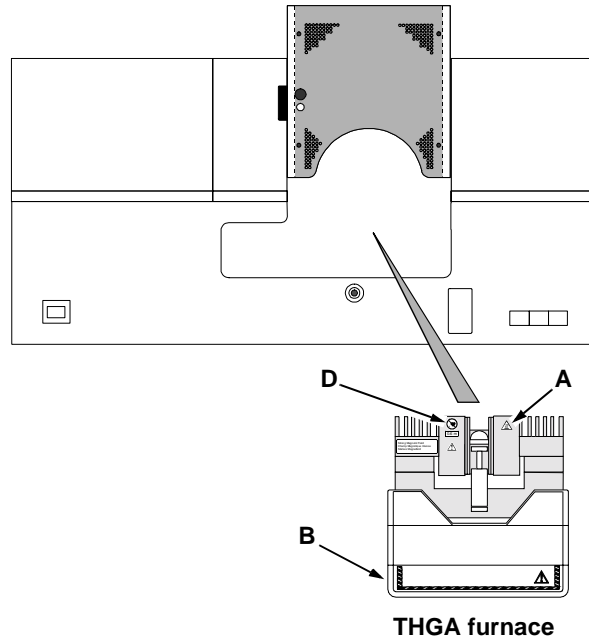
*Anyone wearing a heart pacemaker or having metallic implants should remain at least 0.6 meter away from the furnace – in any direction – while the furnace is operating.*

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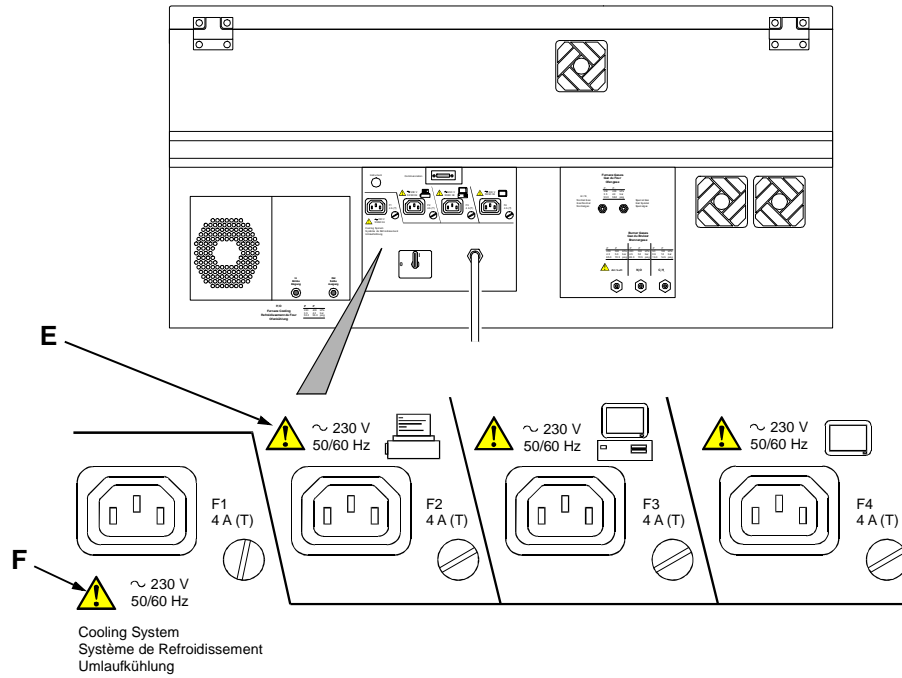
**E**

### **Power outlets**

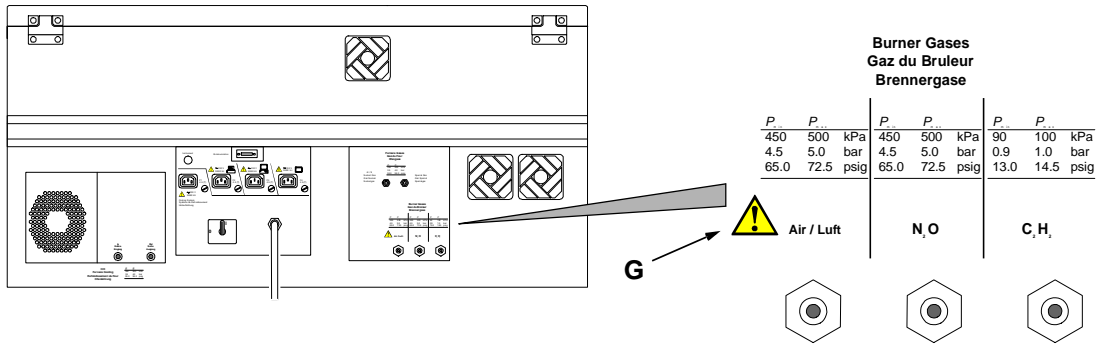
Power outlets for a printer, computer, or monitor.

Use these outlets only when the equipment requires 230 V, 50/60 Hz.

**F**

### **Cooling system outlet**

Connect the cooling system to this power outlet only. Do not connect the cooling system to any other supply. Do not connect other equipment to this power outlet.

G

### Burner Gases

**Never** set the outlet pressure to a value exceeding the maximum pressure ( $P_{max}$ ).

#### Warning: Acetylene – Explosion Hazard

Acetylene can decompose explosively at pressures higher than 103 kPa / 1.03 bar / 15 psig. Always make sure that the acetylene outlet gauge pressure is below this value.

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#### Warning: Flashback Hazard

Perkin Elmer burner systems are designed for use with compressed air. The use of oxygen can cause an explosion in the burner system, and oxygen-enriched air can cause a flashback of the flame.

Never use oxygen or oxygen-enriched air with Perkin Elmer burner systems. Use only compressed air.

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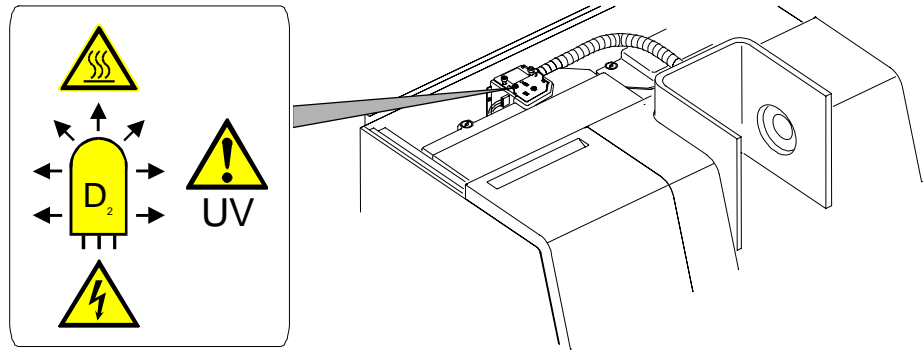
## Warning markings on the deuterium lamp cover

The warning markings below are visible when you open the top cover of the spectrometer.

### **High Temperatures – Risk of Burns**

*The lamp is hot.*

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### **UV Radiation – Risk of Eye Damage**

*The lamp emits intense UV radiation which can damage your eyes.*

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### **Electrical Hazard**

*The deuterium lamp operates at high voltage.*

*To avoid electrical shock, disconnect the line power cord before opening the cover.*

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<b>Contents</b>	<b>page</b>
Introduction . . . . .	2-3
Correct use of analytical instruments . . . . .	2-3
Laboratory hygiene . . . . .	2-4
Laboratory ventilation . . . . .	2-4
Safe handling of chemicals . . . . .	2-5
Waste disposal . . . . .	2-7
Contents of waste containers . . . . .	2-7
Disposing of defective lamps . . . . .	2-7
UV radiation . . . . .	2-8
Magnetic field . . . . .	2-8
Zeeman graphite furnace systems . . . . .	2-8
High temperatures . . . . .	2-9
Burner system . . . . .	2-9
Graphite furnace . . . . .	2-9
Quartz tube atomizer cell . . . . .	2-9
Handling compressed gases . . . . .	2-10
Summary of gas hazards . . . . .	2-10
Identifying cylinders . . . . .	2-10
Storing cylinders . . . . .	2-11
Handling cylinders . . . . .	2-12

<b>Contents</b>	<b>page</b>
Safety practices for flame atomization .....	2-13
Safety interlocks .....	2-13
Safe use of burner gases .....	2-13
Air supply .....	2-14
Nitrous oxide .....	2-14
Acetylene .....	2-15
Safe operation of the flame .....	2-16
The drain system .....	2-17
Flashbacks .....	2-17
Hazards with flame atomization .....	2-18
Safety practices for electrothermal atomization .....	2-19
Safety interlocks .....	2-19
Inert gas .....	2-19
Safety checks .....	2-19
Hazards with electrothermal atomization .....	2-20
Safety practices for FIAS and mercury/hydride systems .....	2-21
FIAS-Cells – To all users who have an MHS-20 in addition to FIAS .....	2-21
Safety checks .....	2-21
Hazards with FIAS and mercury/hydride systems .....	2-22
References for laboratory safety practice .....	2-23

## ***Introduction***

The guides provided with your analytical instrument contain information and warnings that you must follow to ensure safe operation and to maintain the instrument in a safe condition. This advice is intended to supplement, not supersede, the normal safety code of behavior prevailing in the country of operation.

This chapter describes general safety practices for atomic spectroscopy and potential hazards with various atomic absorption techniques. The information provided does not cover every safety procedure that should be practiced. Ultimately, maintenance of a safe laboratory environment is the responsibility of the user and the user's organization.

## ***Correct use of analytical instruments***

Before you install or use your instrument, and in order to get the best results, you should be familiar with all of the instruments in the system and know how to operate them. You should also be aware of the safety procedures in force in your laboratory, especially those concerning atomic spectroscopy instruments. Read the guides supplied with the instruments before you start.

If you use the instrument in a manner not specified in the guides, or if you use it for a purpose other than that intended, you may damage the instrument, or compromise your own, or someone else's, safety.

Analytical instruments should only be operated by persons who are suitably qualified and have received adequate training.

## ***Laboratory hygiene***

- Keep the work area scrupulously clean to avoid contaminating your samples and to maintain a safe working environment. Clean up spilled chemicals immediately and dispose of them properly.
- Do not allow smoking in the work area. Smoking is a source of significant contamination and also a potential route for ingesting harmful chemicals.
- Do not store, handle, or consume food in the work area.

## ***Laboratory ventilation***

Toxic combustion products, metal vapor, ozone, etc., can be generated by the system, depending on the type of analyses being performed.

- You must provide an efficient laboratory ventilation system to remove toxic products generated during instrument operation.

## Safe handling of chemicals

Some chemicals used with the instrument may be hazardous or may become hazardous after completion of an analysis.

The responsible body<sup>1</sup> must take the necessary precautions to ensure that the surrounding workplace is safe and that instrument operators are not exposed to hazardous levels of toxic substances (chemical or biological) as defined in the applicable national, state, and local health and safety regulations and laws. Venting for fumes and disposal of wastes must be in accordance with all national, state, and local health and safety regulations and laws.

- Use, store, and dispose of chemicals in accordance with the manufacturer's recommendations and the applicable national, state, and/or local regulations.
- Wear appropriate eye protection at all times while handling chemicals. Depending on the types of chemicals you are handling, wear safety glasses with side shields, or goggles, or a full-face shield.
- Wear suitable protective clothing, including gloves if necessary, resistant to the chemicals you are handling.
- When preparing chemical solutions, always work in a fume hood that is suitable for the chemicals you are using.
- Perform sample preparation away from the instrument to minimize corrosion and contamination.
- Clean up spills immediately using the appropriate equipment and supplies, such as spill cleanup kits.
- Do not put open containers of solvent near the instrument.
- Store solvents in an approved cabinet (with the appropriate ventilation) away from the instrument.

---

1. Definitions from IEC 1010-1:

**Responsible body:** 'individual or group responsible for the use and maintenance of equipment, and for ensuring that operators are adequately trained.'

**Operator:** 'person operating equipment for its intended purpose.'

**Sodium tetrahydroborate**

Sodium tetrahydroborate is used as a reductant in the mercury cold vapor technique and the hydride-generation technique. Sodium tetrahydroborate solutions are unstable and decompose, releasing hydrogen. Sodium tetrahydroborate also releases hydrogen when it comes into contact with acids.

- Make sure that the work area has an adequate ventilation system to prevent the build-up of explosive hydrogen/air mixtures.
- Use and store sodium tetrahydroborate according to the manufacturer's recommendations.
- Keep sodium tetrahydroborate solutions:
  - Out of direct sunlight.
  - Away from open flames.
  - In an area with an efficient ventilation system.
- Protect sodium tetrahydroborate solutions from temperature variations.
- Handle sodium tetrahydroborate solutions with care since they are corrosive.
- Do not allow smoking in areas where sodium tetrahydroborate is used.
- When you are using sodium tetrahydroborate as the reductant with flow-injection systems, hydrogen is formed when it comes into contact with the acidic carrier solution.
  - Do not loosen or remove any connections when the system is operating.
  - The maximum concentration of sodium tetrahydroborate that you should use with Perkin Elmer flow-injection systems is 0.5% w/v.
  - At the end of the work session, pump deionized water through the sodium tetrahydroborate tubing system to remove all traces of the reductant.

## Waste disposal

### Contents of waste containers

The materials that you collect in waste containers may include small amounts of the substances that were analyzed and other chemicals used in the analyses.

If these materials are toxic, corrosive, or contain organics you may have to treat the collected effluent as hazardous waste. Refer to your local safety regulations for proper disposal procedures.

### Disposing of defective lamps

Hollow cathode lamps and electrodeless discharge lamps contain small quantities of the lamp element in a very pure form. If a lamp containing toxic elements becomes defective or otherwise unusable, you must treat it as hazardous waste and dispose of it accordingly. A licensed company in the chemical waste disposal business can provide lamp disposal in accordance with environmental regulatory requirements. Please note that Perkin Elmer does not take back defective lamps for disposal.

In addition, hollow cathode lamps and deuterium lamps are maintained under reduced pressure. Handle and dispose of them correctly to minimize the implosion risk.

For information, the quantities of analyte material contained in the cathodes of hollow cathode lamps are listed in the table below:

Quantity of material in the cathode	Single-element hollow cathode lamps	Multi-element hollow cathode lamps
less than 5 g	As, Au, B, Ba, Be, Ca, Dy, Er, Eu, Ga, Gd, Ge, Hf, Ho, In, Ir, K, La, Li, Mg, Na, Nd, Pd, Pr, Pt, Re, Rh, Ru, Sc, Se, Sm, Sn, Ta, Tb, Tm, Yb, Y	Na-K, Pt-Ru
5–10 g	Ag, Al, Bi, Cd, Co, Cr, Mn, Mo, Ni, Sb, Si, Sr, Te, V, W, Zn, Zr	Ca-Mg, Ca-Zn, Ag-Au, Sn-Te, Ca-Mg-Zn, Ca-Mg-Al
10–15 g	Cu, Fe, Hg, Nb, P, Pb, Ti, Tl	all other multi-element HCLs

For EDLs, the quantity of analyte material used is much less than the quantities used in HCLs. Typically, only several milligrams of material are placed in the EDL bulb. No Perkin Elmer EDLs (System 1 or 2) contain more than 0.05 g of analyte material.

## ***UV radiation***

You should be aware of the health hazard presented by UV radiation.

- When the instrument is on, do not remove any covers unless specifically instructed to do so in the guide since otherwise you may be exposed to potentially hazardous UV radiation.
- Always wear UV-absorbing eye protection when viewing any of these sources:
  - The flame, especially the nitrous oxide/acetylene flame.
  - The graphite furnace when it is heated to incandescence.
  - Hollow cathode or electrodeless discharge lamps.
  - The deuterium background correction lamp.

## ***Magnetic field***

### ***Zeeman graphite furnace systems***

The electromagnet generates a strong magnetic field inside the furnace during measurement cycles.

There is a weak, accessible, stray magnetic field around the graphite furnace. This stray magnetic field can affect heart pacemakers and metallic implants up to 0.6 meter from the furnace in all directions. This stray field can also affect other instruments, watches, and magnetic storage devices.

- Do not operate the furnace when persons wearing electronic heart pacemakers, or having other metallic implants, are present.
- Anyone wearing a heart pacemaker or having metallic implants should remain at least 0.6 meter away from the furnace – in any direction – while the furnace is operating.

## High temperatures

### **Burner system**

The flame can generate temperatures of up to 2800 °C.

- Do not touch the burner head until it has cooled to room temperature.

### **Graphite furnace**

The THGA graphite furnace can generate temperatures of up to 2600 °C.

- Do not touch any part of the graphite furnace until it has cooled to room temperature.
- Do not attempt to inject a sample into the graphite tube when the furnace is hot. You may be subjected to fumes from the sudden vaporization of the sample, and may damage the pipet.

### **Quartz tube atomizer cell**

The quartz tube atomizer cell (QTA-cell) used for the hydride-generation technique can reach temperatures of up to 1000 °C.

- Do not touch any part of the heating mantle (electrically heated systems), or the cell holder (flame-heated systems), or the QTA-cell until they have cooled to room temperature.

## Handling compressed gases

*Note: The permanent installation of gas supplies is the responsibility of the user and should conform to local safety and building regulations.*

### Summary of gas hazards

Hazards associated with the different gases used in atomic absorption spectrometry are presented in the table below:

Gas	Suffocation	Explosion	Spontaneous Decomposition or Combustion
Air	–	–	–
Argon	x	–	–
Nitrogen	x	–	–
Acetylene (ethyne, C <sub>2</sub> H <sub>2</sub> )	x	x	x
Nitrous Oxide (dinitrogen monoxide, N <sub>2</sub> O)	x	–	x

Contact the gas supplier for a safety data sheet containing detailed information on the potential hazards associated with the gas.

### Identifying cylinders

- Legibly mark cylinders to clearly identify the contents and status (full, empty, etc.). Use the chemical name or commercially accepted name for the gas.

## **Storing cylinders**

- Store cylinders in accordance with the applicable national, state, and/or local regulations and standards.
- When gas cylinders are stored in confined areas, such as a storage room, make sure that ventilation is adequate to prevent toxic or explosive accumulations of gas.  
The storage room should be well ventilated and dry. This is particularly important in confined areas.
- Do not store cylinders near elevators, gangways, or in locations where heavy moving objects may strike or fall against them.
- Use and store cylinders away from exits and exit routes.
- Locate gas cylinders away from heat sources, including heat lamps.  
Compressed gas cylinders should not be subjected to temperatures above 52 °C (125 °F).
- Do not allow ignition sources in the storage area and keep cylinders away from readily ignitable substances such as gasoline or waste, or combustibles in bulk, including oil.
- Store all gas cylinders only in a vertical position, with the valve cap in place, and fastened securely to an immovable bulkhead or a permanent wall.
- If you are storing cylinders outdoors, store them above ground on a suitable floor where they are protected against temperature extremes (including the direct rays of the sun).

### ***Handling cylinders***

- Move cylinders with a suitable hand truck after making sure that the valve cap is securely in place and that the cylinder is properly fastened to the hand truck.
- Use only approved regulators, tubing, and hose connectors. When you connect fittings, keep in mind that left-hand thread fittings are used for fuel gas connections (e.g., acetylene), whereas right-hand thread fittings are used for oxidant and support gas connections (e.g., nitrous oxide, air).
- Arrange gas hoses where they will not be damaged or stepped on and where things will not be dropped on them.
- Do not ‘crack the valve’ or open the valve of an acetylene cylinder before attaching a regulator.
- Do not attempt to refill gas cylinders.
- Check the condition of pipes, hoses, and connectors regularly, and replace any damaged parts.
- Perform periodic gas leak tests at all joints and seals of the gas system by applying an approved gas leak detection solution.

## ***Safety practices for flame atomization***

### ***Safety interlocks***

Perkin Elmer provides a number of safety interlocks on the burner system to monitor gas pressure and check for the proper setup of the burner head, nebulizer, and drain system. In addition a flame sensor checks that the flame is burning.

- Do not attempt to defeat these interlocks; you may compromise your own, or someone else's, safety.

### ***Safe use of burner gases***

The seepage of fuel gas or fuel gas mixture from the burner system, the drain system, the gas control system, or the gas connections constitutes a serious fire hazard.

- Make sure that there are no breaks or leaks in any of these systems and that all the seals are correctly installed and in good condition.
- Regularly check for leaks at joints and seals using an approved leak test solution.
- When you perform any maintenance or installation procedures, follow the instructions in the guide exactly.
- Do not attempt to service the gas control system yourself. A Perkin Elmer service engineer, or similarly authorized and trained person, must perform the work.
- When you shut down the instrument, for example at the end of the working day, shut all the gas lines at the cylinder or regulator valves. Bleed the lines between the regulator and instrument to atmosphere before switching off the ventilation system.

### ***Air supply***

An air compressor is recommended for the air supply to the burner system. Cylinders of compressed air should only be used as an emergency or short-term solution.

- If you use cylinder air, make sure that it is compressed air and not oxygen-enriched air (e.g. breathing air); the use of oxygen-enriched air can cause a flashback of the flame.
- **Never** use oxygen as the oxidant since this can cause an explosion.

### ***Nitrous oxide***

(dinitrogen monoxide, N<sub>2</sub>O)

- Tubing and fittings carrying nitrous oxide must be free of all oil, grease, and organic materials.  
Spontaneous combustion may occur if nitrous oxide comes into contact with these materials.
- Use a double-stage or heated regulator for nitrous oxide to prevent freezing of the diaphragm and a loss of pressure regulation.
- Do not store nitrous oxide in close proximity to flammable gases, such as acetylene or hydrogen.

## Acetylene

- Always use 'Atomic Absorption Grade' acetylene (ethyne) dissolved in acetone (propan-2-one). This is satisfactory for all Perkin Elmer AA spectrometers.
- Do not store acetylene in close proximity to oxidizing gases, such as nitrous oxide.
- Use approved tubing and fittings for acetylene. Never use copper fittings since acetylene forms an explosive compound with copper.
- Prevent acetylene from coming into contact with copper, silver, mercury or gaseous chlorine.
- Periodically check for the presence of acetylene in the laboratory atmosphere, especially near the ceiling.
- For acetylene, set an outlet gauge pressure of between 85 kPa and 100 kPa (0.85–1.0 bar, 12–14.5 psig). Never allow the outlet gauge pressure to exceed 103 kPa (1.03 bar, 15 psig); acetylene can explode spontaneously above this pressure.
- Solvent Carryover.  
If the acetylene cylinder pressure falls to below 600 kPa / 6.0 bar / 87 psig (at 20 °C), some of the solvent used to stabilize the acetylene (such as acetone) may be carried over into the burner. This could influence the characteristics of the burner with a resulting influence on the analytical results.  
Change to a new acetylene cylinder when the cylinder pressure falls to below this value.

### **Safe operation of the flame**

- Before you ignite the flame make sure that:
  - The laboratory fume ventilation system is operating;
  - The burner head is installed correctly;
  - The burner end cap is secured firmly;
  - The fuel and oxidant tubing fittings are properly connected;
  - The burner unit contains the necessary solvent-resistant components if you intend to analyze samples containing organic solvents;
  - The atomizer compartment door is closed;
  - The pH of the liquid in the waste container is greater than pH 10 if you intend to aspirate cyanide solutions.
- **Do not leave the flame unattended.** Always make sure that there is a fire extinguisher near the instrument.
- Never change the gas pressure or shut a gas valve while the flame is burning.
- **Do not allow the burner head slot to become blocked. This can cause a flashback of the flame.**
- Do not place open containers of flammable liquids and solvents near to the flame. Be especially careful with samples that contain highly volatile solvents.
- When you have analyzed samples containing organic solvents, flush all traces of solvent from the burner system.

## ***The drain system***

- Always place the waste container in a well ventilated place underneath the spectrometer, **in full view** while you are working with the spectrometer. This prevents the build-up of potentially hazardous gases, and allows you to see the liquid level.
- Never use a glass waste container.
- **Do not kink, fold, or loop the drain tube.**
- Never place the drain tube directly into a laboratory sink.
- Empty the waste container frequently; especially do not allow the waste container to fill with organic solvent.
- Thoroughly flush the drain system, as described in the guide, when you shut down the system.
- Check the condition of the drain tube regularly, especially if you use organic solvents. Replace the drain tube when it first shows signs of cracking or discoloration.

## ***Flashbacks***

A flashback is an explosion of the premixed gases in the spray chamber. The safety interlocks normally prevent conditions that could cause a flashback, but if a flashback should ever occur:

- **Check that the burner head slot is clean.**
- Make sure that you are using compressed air and not oxygen-enriched air.
- Check the flow spoiler or impact bead for damage, and replace any damaged parts.
- Carefully check and tighten all connections to, and seals on, the burner unit, and check the gas flow settings before you attempt to ignite the flame. Seepage of fuel and oxidant from the spray chamber may cause a flashback.
- Check that the burner head is seated correctly.

## **Hazards with flame atomization**

### **Toxic combustion products:**

Toxic combustion products can be generated by the system, depending on the type of analyses being performed.

- You must provide an efficient laboratory ventilation system to remove toxic products generated during instrument operation.

### **High temperatures:**

The flame can generate temperatures of up to 2800 °C.

- To avoid serious burns, never touch the burner head until it has cooled to room temperature.

### **UV radiation:**

The flame, especially the nitrous oxide/acetylene flame, emits ultraviolet radiation.

- Keep the atomizer compartment door closed when the flame is burning and never directly view the flame unless you are wearing UV-absorbing glasses.

### **Acetylides:**

If you have aspirated high concentrations of copper, silver, or mercury solutions into an acetylene flame, unstable acetylides may have formed in the spray chamber. If permitted to dry, these compounds may explode.

- Aspirate solution continuously to prevent residues from drying.
- Thoroughly flush the spray chamber and drain system with water immediately after such an analysis.

### **Cyanide solutions:**

If you intend to aspirate cyanide solutions, make sure that the pH of the liquid in the waste container is greater than pH 10; toxic hydrogen cyanide gas is formed when cyanides contact acidic solutions.

- Never allow solutions containing cyanides to mix with acidic solutions.

## Safety practices for electrothermal atomization

### Safety interlocks

Perkin Elmer provides a number of safety interlocks on the graphite furnace (electrothermal atomization) system that monitor the gas pressure and the temperature of certain system components.

- Do not attempt to defeat these interlocks; you may compromise your own, or someone else's, safety.

### Inert gas

The graphite furnace requires a supply of inert gas. Perkin Elmer recommends **argon**.

Although nitrogen can be used, it is not recommended for the inert gas supply. At furnace temperatures above 2300 °C, nitrogen reacts with graphite to form toxic cyanogen gas (CN)<sub>2</sub>.

### Safety checks

- Before you start an analysis, make sure that:
  - The laboratory fume ventilation system is operating;
  - The inert gas supply is connected and set to the correct pressure;
  - The cooling water supply is turned on;
  - The autosampler waste bottle is empty.

## **Hazards with electrothermal atomization**

### **Toxic products:**

Toxic metal vapors, ozone, etc., can be generated by the system, depending on the type of analyses being performed.

- You must provide an efficient laboratory ventilation system to remove toxic products generated during instrument operation.
- If you use nitrogen as the inert gas, toxic cyanogen gas (CN)<sub>2</sub> will be produced when the furnace temperature is above 2300 °C.

### **High temperatures:**

The THGA graphite furnace can generate temperatures of up to 2600 °C.

- To avoid serious burns, never touch the graphite components until the furnace has cooled to room temperature.
- Do not attempt to inject a sample into the graphite tube when the furnace is hot. You may be subjected to fumes from the sudden vaporization of the sample, and may damage the pipet.

### **UV radiation:**

The graphite furnace emits ultraviolet radiation when it is heated to incandescence.

- Never directly view the graphite furnace unless you are wearing UV-absorbing glasses.

### **Strong magnetic field:**

The electromagnet generates a strong magnetic field inside the THGA furnace during measurement cycles.

- Do not operate the furnace when persons wearing electronic heart pacemakers, or having other metallic implants, are present.
- Anyone wearing a heart pacemaker or having metallic implants should remain at least 0.6 meter away from the furnace – in any direction – while the furnace is operating.

## Safety practices for FIAS and mercury/hydride systems

### **FIAS-Cells – To all users who have an MHS-20 in addition to FIAS**

- Never use the FIAS-cell for the MHS-20.
- The smaller diameter of the FIAS-cell can cause an increase in back pressure. This could force the acidic reaction mixture back into the sodium tetrahydroborate storage container where large quantities of hydrogen would be produced.

### **Safety checks**

- Before you start an analysis, make sure that:
  - The laboratory fume ventilation system is operating;
  - The tubing in the fluid system is not constricted since this could lead to a pressure build-up in the system;
  - The drain tubing is inserted in the waste container and that this container is empty.
- If you are using a flame to heat the QTA-cell or if you intend to perform flame-FIAS analyses, refer to the information under *Safety Practices for Flame Atomization* before you ignite the flame.

## **Hazards with FIAS and mercury/hydride systems**

### **Toxic products:**

Toxic metal vapors, etc., can be generated by the system, depending on the type of analyses being performed.

- You must provide an efficient laboratory ventilation system to remove toxic products generated during instrument operation.

### **High temperatures:**

The QTA-cell used for the hydride-generation technique can reach temperatures of up to 1000 °C.

- Do not touch any part of the heating mantle (electrically heated systems), or the cell holder (flame-heated systems), or the QTA-cell until they have cooled to room temperature.

### **Hazardous chemicals:**

Some of the chemicals required for analyses are corrosive and/or toxic.

- Refer to the information under *Safe Handling of Chemicals* and to the specific warnings in the documentation provided with the system before you start an analysis.

## References for laboratory safety practice

- Bretherik, L., *Bretherik's Handbook of Reactive Chemical Hazards*, 4th ed., Butterworth & Co. Ltd., London, UK, 1990.
- Bretherik, L., *Hazards in the Chemical Laboratory*, 3rd ed., Royal Society of Chemistry, London, UK, 1981.
- *CRC Handbook of Laboratory Safety*, 3rd ed., Furr, K., ed., The Chemical Rubber Co. Press, Florida, USA, 1990.
- *Prudent Practices for Handling Hazardous Chemicals in Laboratories*, National Research Council, National Academic Press, Washington D.C., USA, 1981.
- Sax, N., ed., *Dangerous Properties of Industrial Materials*, 7th ed., Van Nostrand Reinhold, New York, 1989.
- Compressed Gas Association, Inc., Arlington, VA 22202, USA, various publications.
- Data Sheets provided by chemical manufacturers, e.g.:
  - USA, Material Safety Data Sheets (MSDS),
  - FRG, DIN-Sicherheitsblätter,
  - UK, Hazard Data Sheets.

<b>Contents</b>	<b>page</b>
Before you install the spectrometer .....	3-3
Moving and reinstalling the spectrometer .....	3-4
Preparing the spectrometer to be moved .....	3-4
Moving the spectrometer .....	3-5
Reinstalling the spectrometer .....	3-5
Connecting the burner gases .....	3-6
Connecting the furnace gases .....	3-9
Connecting the drain system to the burner .....	3-14
Installing the furnace autosampler .....	3-14
Installing the cooling system .....	3-15
Electrical connections .....	3-18

## **Before you install the spectrometer**

### **When the spectrometer arrives**

- Check the exterior of the shipping crate for obvious signs of damage. If the crate has been damaged, file a claim with the authorized carrier immediately and inform your nearest Perkin Elmer office.
- Contact your local Perkin Elmer office to arrange for the installation.

### **Before the service engineer arrives**

- You must prepare a suitable site for the spectrometer and any accessories. See '*Laboratory Requirements*', and the guides for any accessories.
- You must make sure that the fume ventilation system and the necessary gas supply lines and regulators are installed.
- Although you should not perform the initial installation yourself, we recommend that you read the installation instructions to learn about the spectrometer and its requirements before you start using it. These instructions are also provided in case you need to alter any of the connections or move the instrument.

### **When the service engineer arrives**

- The service engineer will remove the spectrometer from the crate and check that all parts have been delivered. If anything is missing or damaged, the service engineer will make sure that the respective parts are supplied.
- The service engineer will remove the shipping locks, perform pre-installation checks, install the spectrometer, and make sure that everything functions correctly.

## **Moving and reinstalling the spectrometer**

We recommend that only a Perkin Elmer service engineer or a Perkin Elmer trained and authorized person moves and reinstalls the spectrometer. Should you attempt to do this yourself, Perkin Elmer cannot be held responsible for any damage or malfunctioning of the spectrometer. The instructions here are primarily intended for occasions where you may want to move the spectrometer **short** distances to a new location within the laboratory. Read the instructions carefully and be sure you understand them thoroughly before you attempt to move the spectrometer. Read the detailed information given in the relevant sections of this chapter before you reinstall the spectrometer. If you intend to transport the spectrometer over a longer distance, contact your local Perkin Elmer office since internal shipping locks must be set via the service software, inside the spectrometer.

### ***Preparing the spectrometer to be moved***

1. Close the AA WinLab application, then shut down the Windows operating system.
2. Switch off the spectrometer, computer and all the accessories.
3. Disconnect all the electrical cables.
4. Remove the furnace autosampler from the spectrometer (*see 'Removing the autosampler from the spectrometer' in the 'Graphite Furnace User's Guide'*).
5. Disconnect the burner gas supplies.
6. Disconnect the drain system from the burner.
7. Disconnect the furnace gas supplies.
8. Prepare the computer and other system components for transport as described in their own guides.
9. Attach the original handles securely. Only use the original handles and screws.

*Make sure that the handles are properly secure before you move the spectrometer.*

## Moving the spectrometer

**Warning: Risk of injury through lifting**

W1.4

The instrument weighs approximately 190 kg.

Improper lifting can cause painful and sometimes permanent back injury.

- Use proper lifting techniques and always use the handles to lift or move the instrument.

**Caution: Risk of damage to the instrument**

The spectrometer is a precision optical instrument. Move the spectrometer very carefully and protect it from vibrations and jolts. Do not tilt the spectrometer.

If the spectrometer must be tilted or mechanically lifted, contact your Perkin Elmer service representative before moving the spectrometer. The shipping locks must be properly secured to prevent any damage occurring if the spectrometer is tilted or jolted.

Carefully lift the spectrometer onto a movable table to transport it to the new location.

## Reinstalling the spectrometer

Reinstall the spectrometer by following exactly the procedures described in this chapter, as summarized in the steps below.

1. Prepare the new site for the spectrometer; see 'Laboratory Requirements'.
2. Remove the handles. Save the handles and screws.  
Insert the plastic plugs provided (B015-3803) into the screw holes.
3. Connect the burner and furnace gas supplies.
4. Connect the drain system to the burner.
5. Install the furnace autosampler.
6. Connect the cooling system for the graphite furnace.
7. Make sure that the main power switches of all the instruments are in the off position, then connect the electrical cables.

## Connecting the burner gases

The connectors for the burner gases are located on the rear panel of the spectrometer.

### **What you need to provide:**

- Suitable supplies of acetylene, air, and nitrous oxide. The gases must conform with the requirements listed in 'Laboratory Requirements'.
- Suitable pressure regulators and on/off valves near to the spectrometer for the gas supplies.

The gas supply systems must comply with your local safety regulations. Refer also to the information given under 'Safety Information'.

### **Warning: Flashback Hazard**

W5.9

*Perkin Elmer burner systems are designed for use with compressed air. The use of oxygen can cause an explosion in the burner system, and oxygen-enriched air can cause a flashback of the flame.*

- *Never use oxygen or oxygen-enriched air with Perkin Elmer burner systems. Use only compressed air.*

### **What is provided with the spectrometer:**

- Gas hose assemblies for acetylene, air, and nitrous oxide.  
The fittings on the gas hose assemblies are suitable for connection to most commercially available pressure regulators. These are R 3/8" standard threaded fittings. The fuel (acetylene) fitting has a left-hand thread, while the oxidant (air, nitrous oxide) fittings have right-hand threads.  
If you require an adapter for your regulator, contact the gas supplier or the regulator manufacturer, or your local Perkin Elmer office.

**To connect the gases**

**Air**

1. Make sure that the air supply is turned off.
2. Use the black hose assembly to connect a source of clean, dry, and filtered air to the fitting labelled **Air/Luft** on the rear panel of the spectrometer.

**Nitrous oxide**

1. Make sure that the nitrous oxide supply is turned off.
2. Use the blue N<sub>2</sub>O hose assembly to connect a source of clean, dry nitrous oxide to the fitting labelled **N<sub>2</sub>O** on the rear panel of the spectrometer.

**Acetylene**

1. Make sure that the acetylene supply is turned off.
2. Use the red C<sub>2</sub>H<sub>2</sub> hose assembly to connect a source of clean, dry acetylene to the fitting labelled **C<sub>2</sub>H<sub>2</sub>** on the rear panel of the spectrometer.

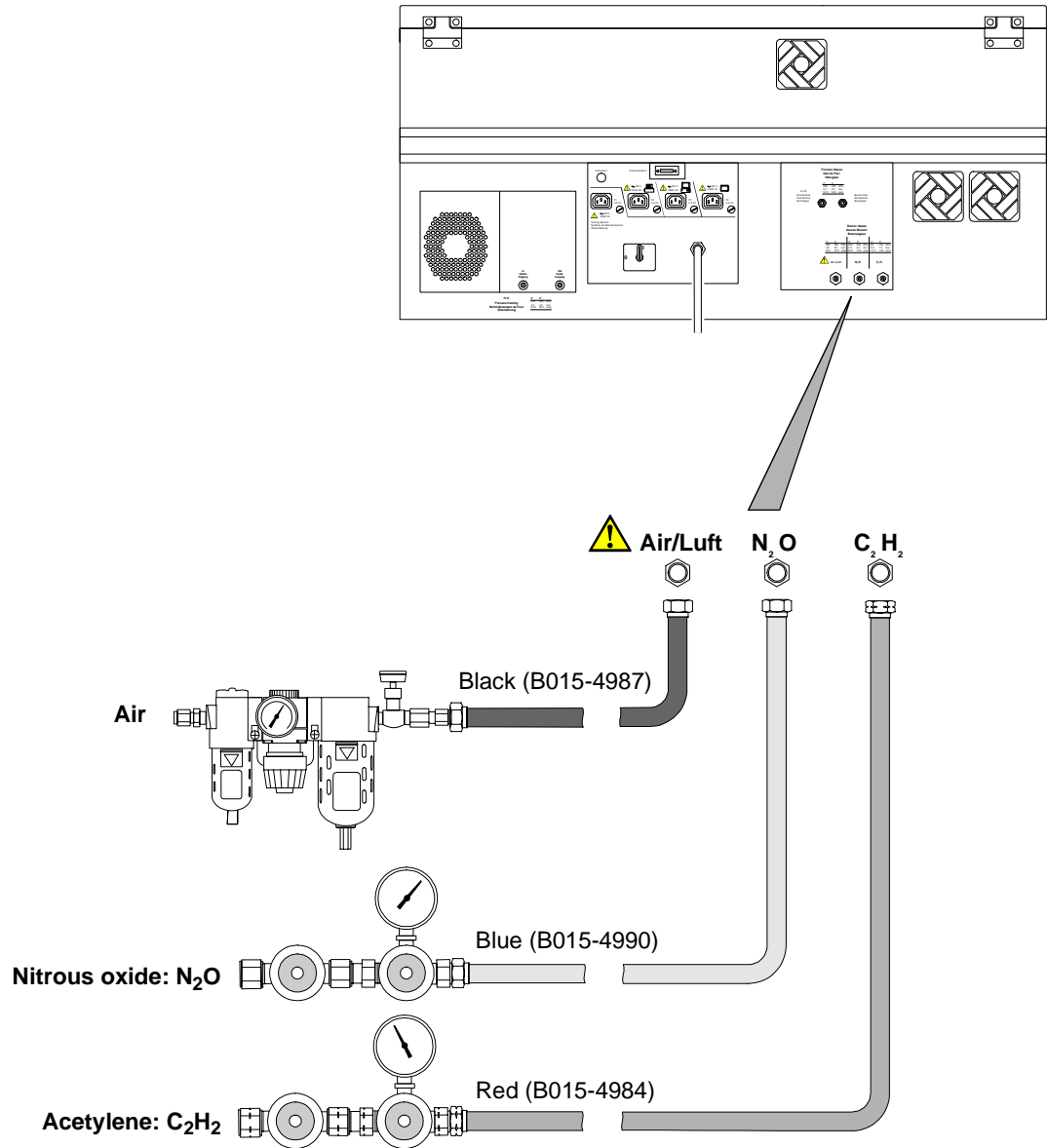
Figure 3-1. Connecting the Burner Gases

## Connecting the furnace gases

The connectors for the furnace gases are located on the rear panel of the spectrometer.

**Note:** *The furnace will operate correctly only if there is gas pressure at both the Normal Gas and Special Gas inlets. If you are using only an inert gas you must connect the gas to both inlets using the Y-connector.*

### **What you need to provide:**

- A suitable supply of argon (or nitrogen) and if necessary a suitable supply of a special gas. The gases must conform with the requirements listed in *Laboratory Requirements*.
- Suitable pressure regulators and on/off valves near to the spectrometer for the inert and special gas supplies.
- A tube cutter for the special gas tube. This is an optional accessory; Part No. 0992-3183.

The gas supply systems must comply with your local safety regulations. Refer also to the information given under '*Safety Information*'.

### **What is provided with the spectrometer:**

- About 10 m of special gas tube, B050-8264.
- A Y-connector with automatic push-fit connectors, 0992-0179.
- Two screw connectors, R 1/4" for the gas regulator. These are not supplied with all instrument versions.

**To connect argon (or nitrogen) only**

Refer to Figure 3-2 on the following page.

1. Make sure that the argon (or nitrogen) supply is turned off.
2. Use the special tube cutter to cut two, 10 cm lengths of tube from the special gas tube. The ends of the tubes must be cut squarely.
3. Connect the long gas tube to the argon (or nitrogen) pressure regulator. Use the R ¼" screw connector provided or use a connector or adapter suitable for the gas regulator. Make sure that all the connections are secure.
4. Pass the gas tube to the rear of the spectrometer. Make sure that it cannot be damaged or strained.
5. Push the two short tubes and the long tube into the connectors on the Y-connector as shown on the following page.
6. Push the free ends of the two short gas tubes into the **Normal Gas** and **Special Gas** connectors on the manifold as shown.  
Push the tubes into the connectors until the tubes butt up to the end of the connector.

**Note:** *To disconnect the gas tube from a connector:  
Carefully push the black locking ring of the connector into the connector and pull the tube out of the connector.*

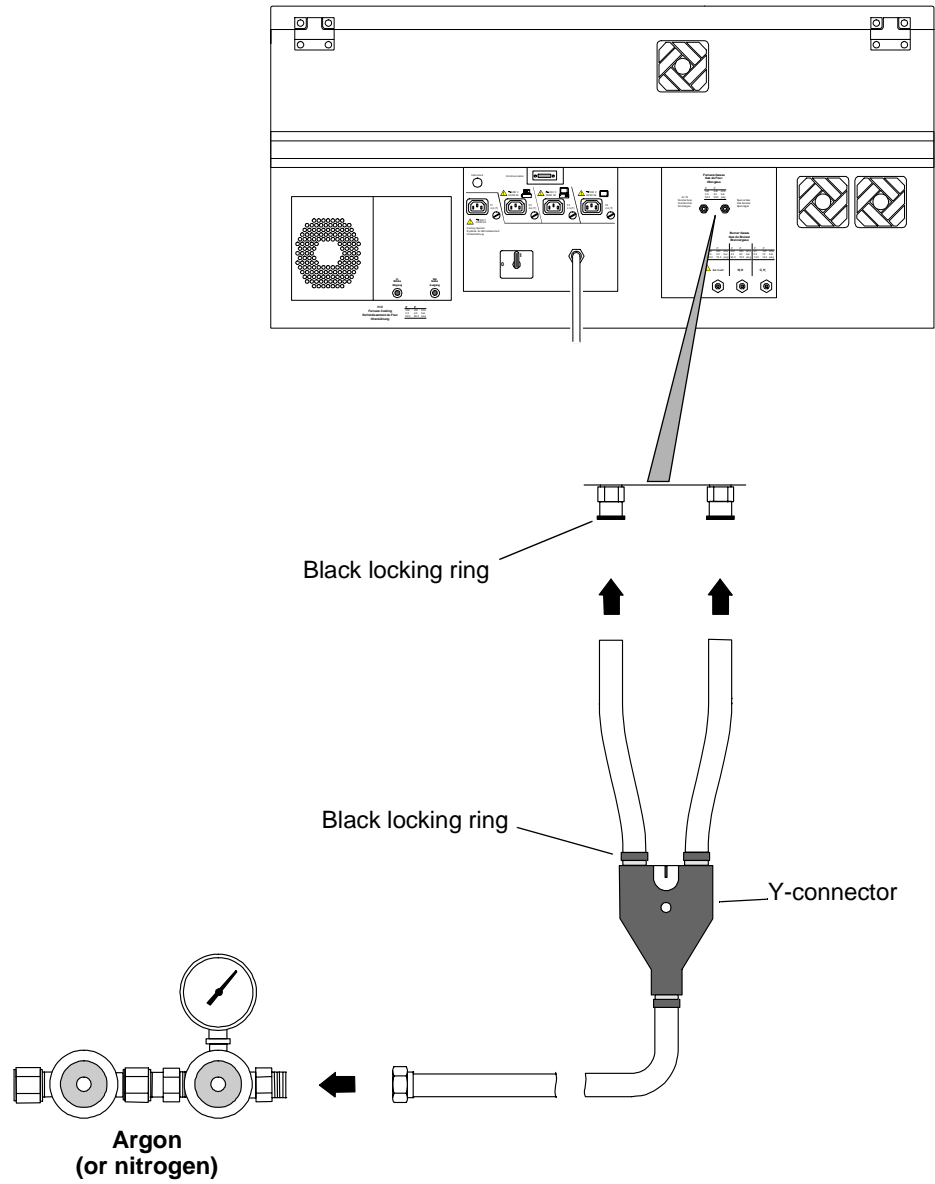
Figure 3-2. Connecting Argon (or Nitrogen) only

**To connect both argon and a special gas**

Refer to Figure 3-3 on the following page.

1. Make sure that the gas supplies are turned off.
2. Use the special tube cutter to cut the special gas tube into two pieces for the two gas supplies. The ends of the tubes must be cut squarely.
3. Connect the gas tubes to the argon (or nitrogen) and special gas pressure regulators. Use the R 1/4" screw connectors provided or use a connector or adapter suitable for the gas regulator. Make sure that all the connections are secure.
4. Pass the gas tubes to the rear of the spectrometer. Make sure that they cannot be damaged or strained.
5. Push the free ends of the gas supply tubes into the **Normal Gas** and **Special Gas** connectors on the manifold as shown.  
Push the tubes into the connectors until the tubes butt up to the end of the connector.

***Note:** To disconnect the gas tube from a connector:*

*Carefully push the black locking ring of the connector into the connector and pull the tube out of the connector.*

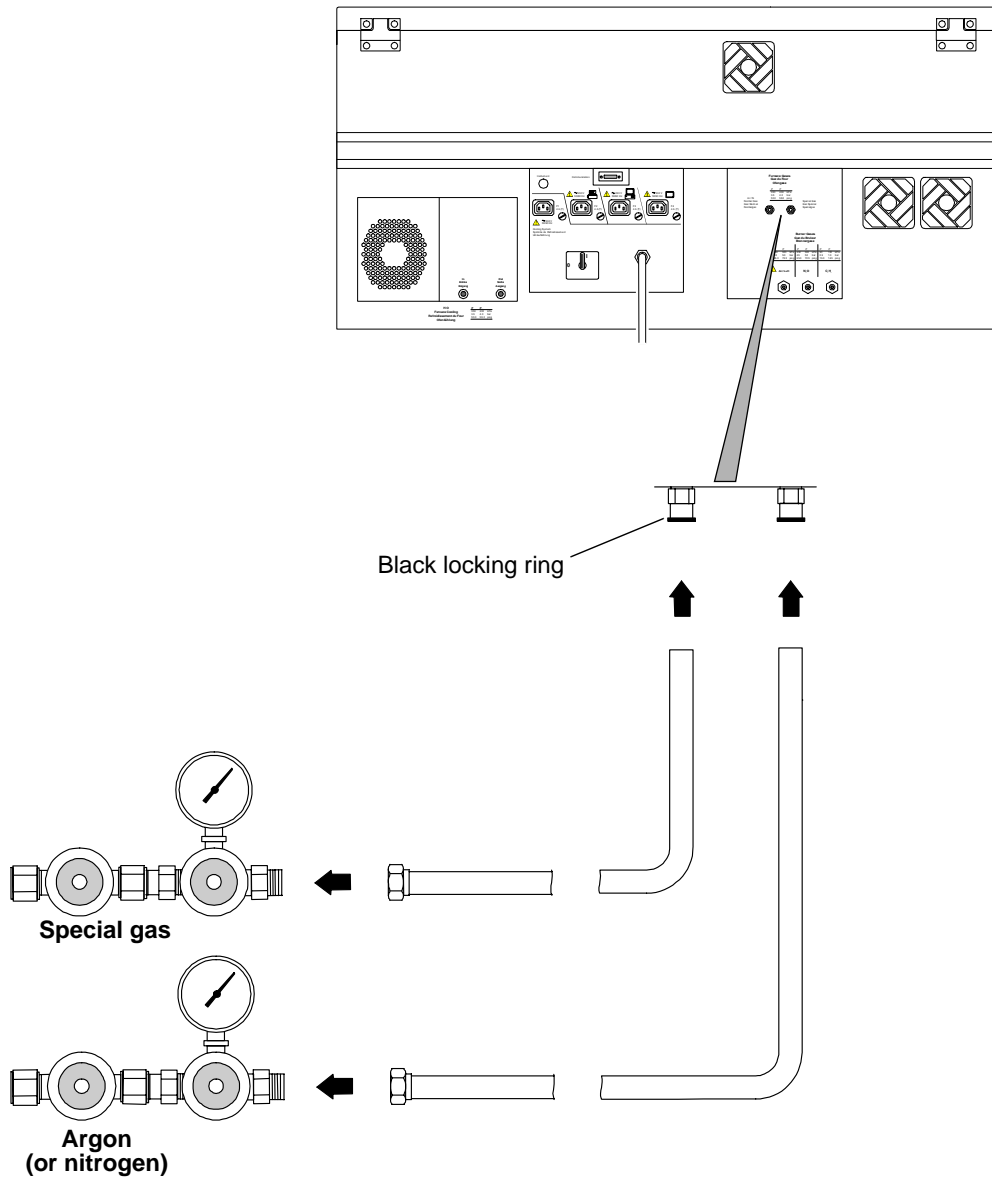
Figure 3-3. Connecting Argon (or Nitrogen) and a Special Gas

### **Connecting the drain system to the burner**

Procedures for connecting the drain system to the burner are described in '*Maintaining the drain system*' in the '*Burner System User's Guide*'.

### **Installing the furnace autosampler**

Procedures for demounting and reinstalling the furnace autosampler are described in '*Installing the furnace autosampler*' in the '*Graphite Furnace User's Guide*'.

## Installing the cooling system

A recirculating cooling system is provided with the spectrometer system. This system includes a pressure-sensitive safety switch that switches off the pump if the internal pressure in the circulatory system increases to above a safe value (see 'Cooling the furnace' in the 'Graphite Furnace User's Guide').

---

**Caution: Risk of damage to the instrument from high water pressure**

High water pressure may cause the coolant tubes on the furnace to burst. The cooling system includes a pressure-sensitive safety switch that switches off the pump if the pressure becomes too high for any reason. The pressure in the main laboratory water supply may be too high for safety.

- Never connect the spectrometer system directly to the main laboratory water supply. Always use the cooling system.
- 

## Connecting and disconnecting the water hoses

The inlet and outlet hoses are permanently attached to the cooling system. They have automatic connectors which are marked with the direction of the water flow.

1. Place the cooling system in a convenient position **on the floor** underneath the spectrometer. Do not place the cooling system on the same bench as the spectrometer since vibrations can adversely influence the spectrometer.

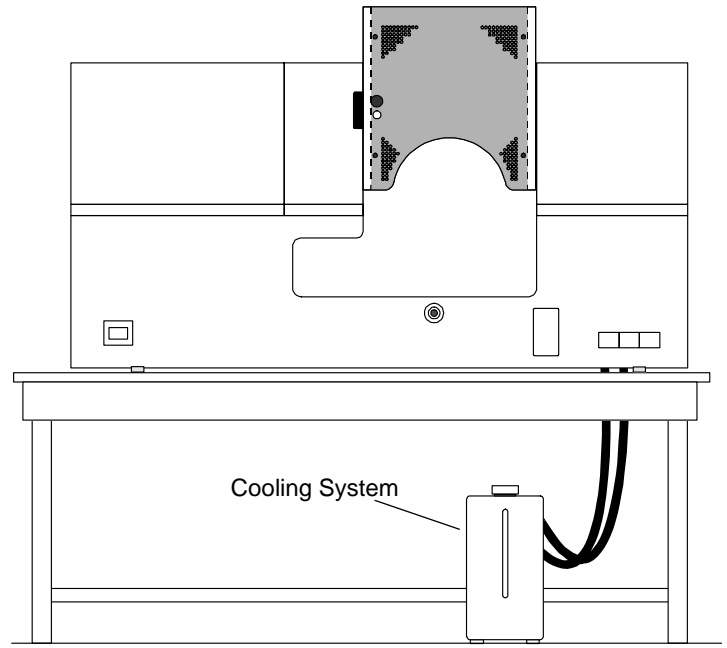
Make sure that:

- You can reach the on/off switch at the rear.
  - There is space at the rear and sides of the cooling system for air to circulate freely.
  - There is space to remove the stopper and fill the system.
  - You can see the water level indicator.
2. Pass the two hoses from the cooling system to the rear of the spectrometer. Make sure they are not strained.
  3. Push the hose connectors onto the correct fittings on the spectrometer as shown opposite.  
The connectors are spring loaded. Push them straight onto the fittings until they click into place.


**Note:** To disconnect the water hoses:

Carefully pull the knurled locking ring down slightly, to release the lock, and pull the connector off the fitting on the spectrometer.

Locating the cooling system

Connecting the hoses

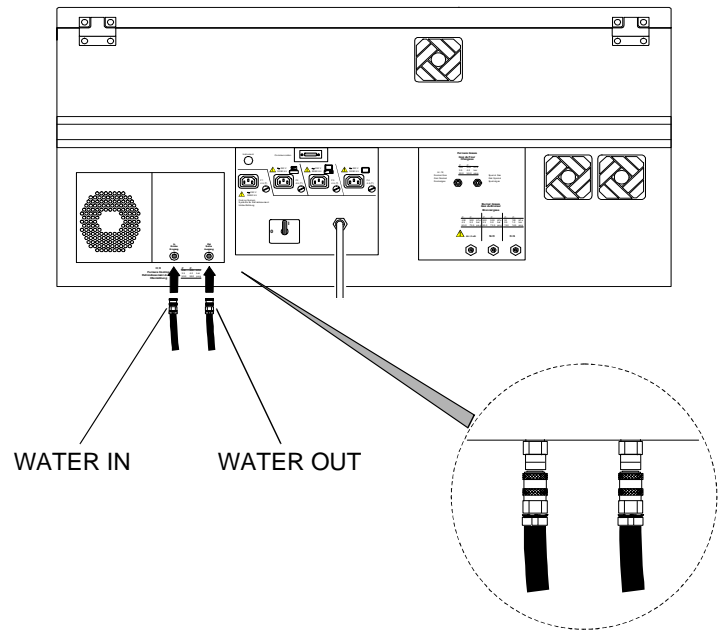
### **Filling the cooling system**

The cooling system is self-priming; air in the system is removed automatically as the cooling system operates. The coolant is a 1:10 solution of glycerine in water (1 part glycerine plus 9 parts water).

**Note:** The cooling system contains 0.5 L glycerine when it leaves the factory. (See 'Cooling system maintenance procedures' in the 'Graphite Furnace User's Guide' for the procedure to follow when you empty and refill the cooling system. Do not use antifreeze.)

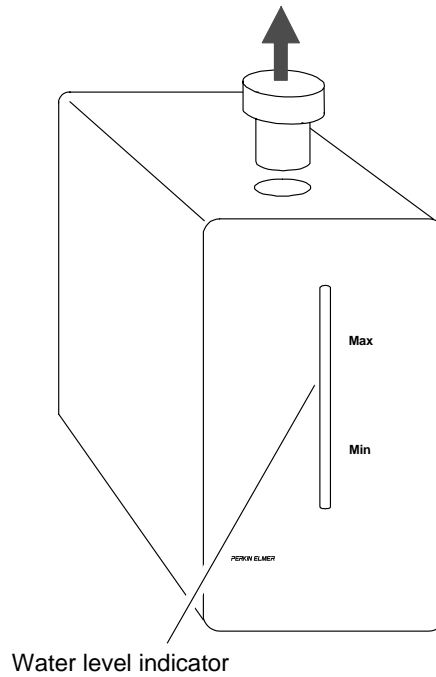1. Remove the stopper from the top of the cooling system – twist and pull.
2. Fill the system to the **Max** mark with deionized water or sediment free drinking water – about 4.5 L.
3. Replace the stopper securely.

## **Electrical connections**

**Warning: Electrical Hazard**

W2.3

- To prevent potential injury to yourself and damage to the instrument, switch OFF all instruments in the system and disconnect them from the line power supply before you perform the procedures described below.

**Read the following instructions very carefully before connecting any cables.**

See 'Laboratory Requirements' for the specification of the line power supply.

### **Summary**

1. Make sure that the spectrometer and the cooling system are switched off.
2. Switch the circuit breaker on the rear panel of the spectrometer to the off position (marked ○).
3. Switch off the computer and all other instruments that you will connect to the spectrometer or computer. See the guides for the individual instruments for the procedure to use.
4. Disconnect the line power cords of the spectrometer and all other instruments from the line power supply.
5. Connect the instrument control cables and power cords as shown on the following page.
6. When you have completed all connections, switch on the system (see 'Switching on the system', page 4-3).

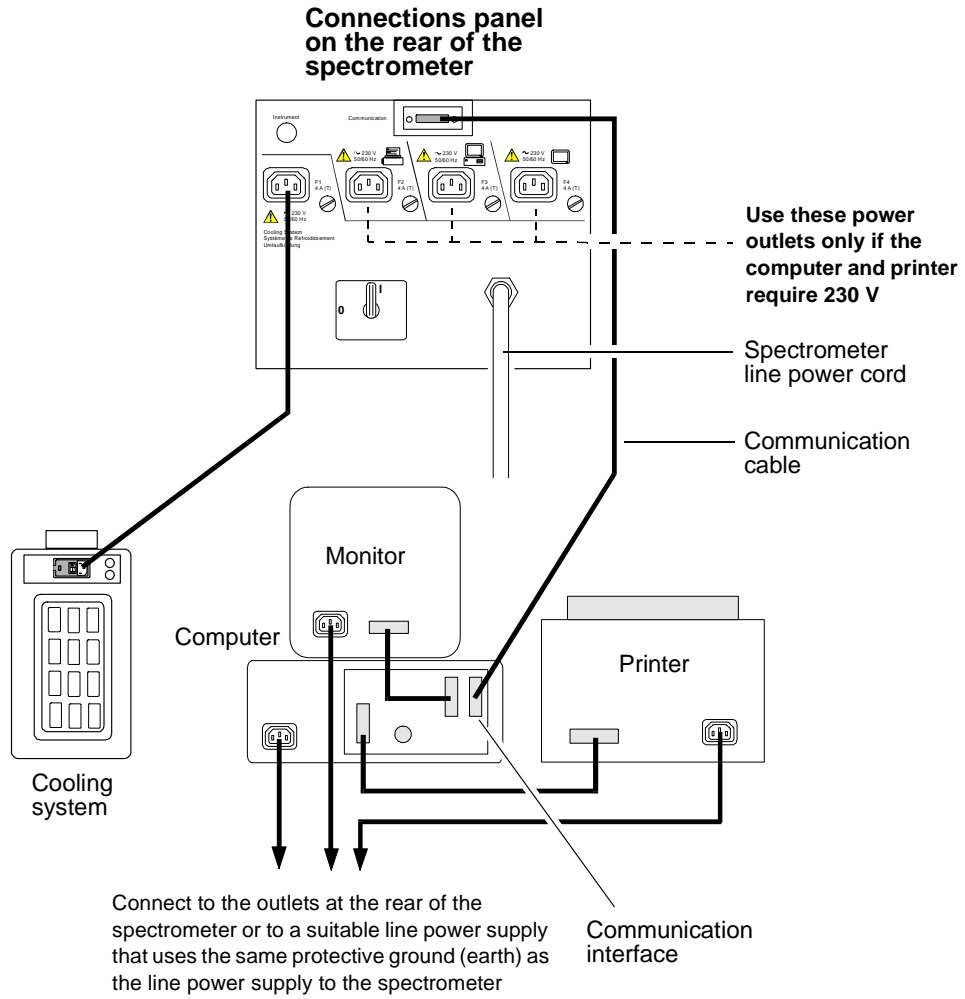
Figure 3-4. Electrical Connections

# Preparing the System for Analyses

---

# 4

<b>Contents</b>	<b>page</b>
Switching on the system .....	4-3
Installing lamps .....	4-7
Installing sample trays .....	4-11
Shutting down the system .....	4-12
Emergency shutdown .....	4-14

## Switching on the system

1. Read the safety information at the front of this guide before you operate the system.
2. Make sure that the spectrometer system and other ancillary instruments are correctly installed and connected to their power supplies.  
Make sure that the coolant hoses are connected correctly to the spectrometer. Where applicable, make sure that the cooling system is filled to the **Max** mark (see 'Filling the cooling system', page 3-17).
3. Switch on the fume ventilation system for the work area.
4. Turn on the burner gases and adjust the outlet gauge pressures to the recommended values (see Table 4-1 on page 4-5).
5. Turn on the inert (and special) gas for the furnace and adjust the outlet gauge pressures to the recommended values (see Table 4-2 on page 4-6).

**Note:** You must always turn on the air supply since air pressure is required for the internal pneumatic system. If you intend to work with only one atomizer, you only need to additionally turn on the gas supplies for that atomizer.

6. Turn on the cooling water to the system:  
Switch on the circulatory cooling system by pressing the power switch located on the rear panel.  
*You can leave this switch permanently in the on position since power is supplied by the spectrometer.*
7. Switch the circuit breaker on the rear panel of the spectrometer to the on position (marked I).  
*You can leave this switch permanently in the on position during routine operation; you switch the spectrometer system on and off via the operational on/off switch located on the front panel.*
8. Switch on the computer.  
If your computer has a power management system, switch this off.
9. Install a Lumina hollow cathode lamp in any holder in the lamp compartment (procedure: page 4-7).  
*A hollow cathode lamp must be installed in at least one holder before you switch on the spectrometer.*
10. Switch on the spectrometer system with the operational on/off switch located on the front panel.

11. If you need the diluter, autosampler, or flow injection system, switch on the respective systems.
12. If you have changed the configuration of the system, such as installed a different autosampler or added or removed any accessories, use the **Reconfigure** utility to reconfigure the software (*see the Online Help in AA WinLab*).
13. Start AA WinLab: In the AA WinLab group, double click on the AA WinLab icon.
14. Set up the various parts of the system:
  1. Install any further lamps that you require for the analyses (*procedure: page 4-7*).
  2. Set up the burner system (*see 'Setting up the system for analyses using the flame technique' in the 'Burner System User's Guide'*).
  3. Set up the graphite furnace system (*see 'Setting up the system for analyses using the furnace technique' in the 'Graphite Furnace User's Guide'*).

Procedures for shutting down the system are described on page 4-12.

In an emergency, you can shut down the system as described on page 4-14.

Table 4-1. Required Pressures for the Burner Gases

Gas	Outlet Gauge Pressure					
	kPa		bar		psig	
	$P_{min}$	$P_{max}$	$P_{min}$	$P_{max}$	$P_{min}$	$P_{max}$
Acetylene (C <sub>2</sub> H <sub>2</sub> )	90	100	0.9	1.0	13.0	14.5
Air	450	500	4.5	5.0	65.0	72.5
Nitrous oxide (N <sub>2</sub> O)	450	500	4.5	5.0	65.0	72.5

- Set the outlet gauge pressure for each gas to a value just below the maximum value ( $P_{max}$ ).
- Always set the same outlet gauge pressure for air and nitrous oxide.
- **Never** set the outlet gauge pressure to a value exceeding the maximum pressure ( $P_{max}$ ).
- Always turn on the air supply since air pressure is required for the spectrometer's pneumatic system.

**Warning: Acetylene – Explosion Hazard**

W5.7

Acetylene can decompose explosively at pressures higher than 103 kPa/ 1.03 bar/15 psig.

- Always make sure that the acetylene outlet gauge pressure is below this value.

Table 4-2. Required Pressures for the Furnace Gases

Gas	Outlet Gauge Pressure					
	kPa		bar		psig	
	$P_{\min}$	$P_{\max}$	$P_{\min}$	$P_{\max}$	$P_{\min}$	$P_{\max}$
Argon	350	400	3.5	4.0	50.0	58.0
Nitrogen	350	400	3.5	4.0	50.0	58.0
Special	350	400	3.5	4.0	50.0	58.0

- Set the outlet gauge pressure for each gas to a value just below the maximum value ( $P_{\max}$ ).
- **Never** set the outlet gauge pressure to a value exceeding the maximum pressure ( $P_{\max}$ ).

**Warning: Toxic Gases**

W6.1

If you use nitrogen as the inert gas, toxic cyanogen gas (CN)<sub>2</sub> will be produced when the furnace temperature is above 2300 °C.

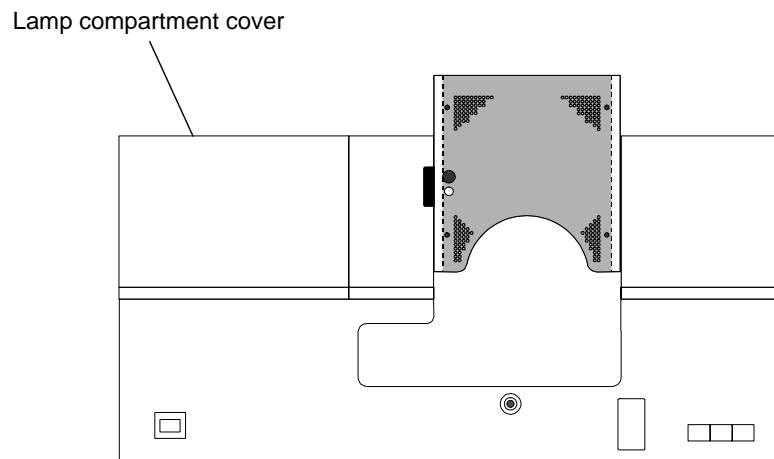
- Make sure that the main laboratory fume ventilation system removes all gases and fumes emitted by the furnace.

## Installing lamps

### Lamp compartment

The lamp compartment is located on the left-hand side of the spectrometer, beneath a hinged cover. The compartment is easily accessible from the front when the cover is open.

The spectrometer is designed to operate with Perkin Elmer Lumina hollow cathode lamps (HCLs) and coded electrodeless discharge lamps (EDLs).

The lamp compartment has holders for up to eight lamps.

You can install Lumina HCLs in all eight holders. The holders incorporate the receptacles for connection of the Lumina lamps.

You can install EDLs in holders 1–4 only. The sockets for the power supplies to the EDLs are located on the left-hand side of the lamp compartment.


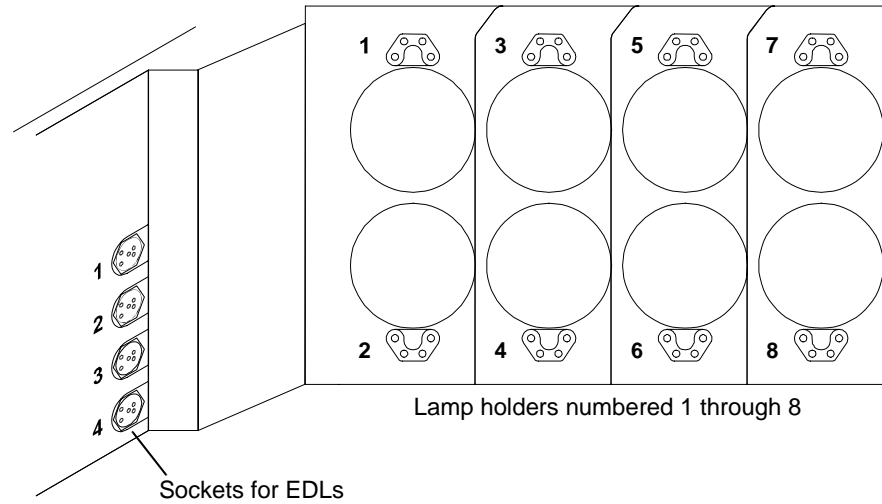
Figure 4-1. Lamp Holders in the Lamp Compartment

### Installing and connecting lamps
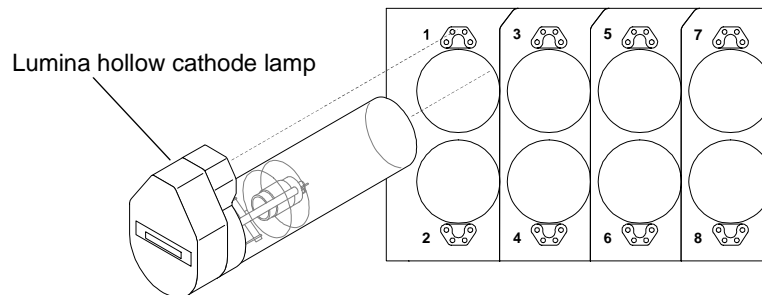
#### HCLs

**Note:** A hollow cathode lamp must be installed in at least one holder before you switch on the spectrometer.


1. Open the lamp compartment cover by lifting it at the front and swinging it back.
2. Slide the Lumina lamp for the required element into the selected lamp holder and make sure that the integrated socket slides fully home over the receptacle on the lamp holder.
3. Close the lamp compartment cover.


**EDLs**

You must use the lamp holders numbered 1 through 4 for EDLs.

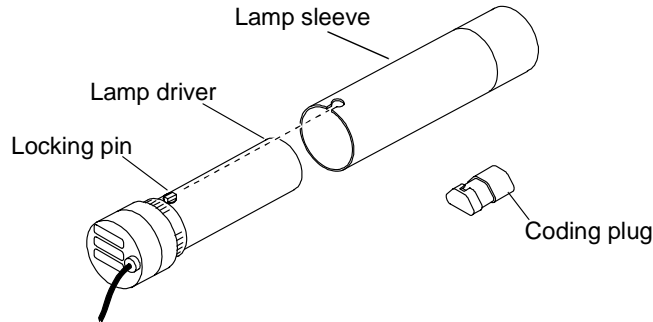1. Open the lamp compartment cover by lifting it at the front and swinging it back.
2. If necessary, assemble the lamp by sliding the lamp driver into the lamp sleeve for the required element. Depress the locking pin as the sleeve slides into place and make sure that the locking pin extends through the locating hole in the sleeve.

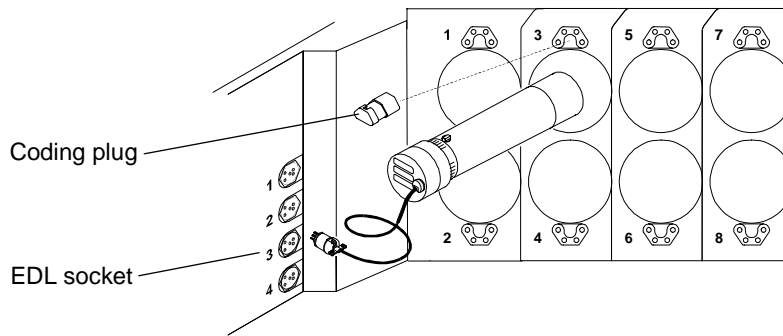3. Slide the lamp carefully into the selected lamp holder until it butts up against the end-stop inside the holder.
4. Insert the coding plug over the receptacle on the lamp holder.
5. Insert the lamp connector into the correct EDL socket at the left-hand side of the lamp compartment. Use the receptacle that has the same number as the lamp holder.
6. Close the lamp compartment cover.

### **Removing lamps**

Do not unplug and remove lamps while they are switched on.

1. On the **Toolbar**, click on **Lamps**. The Align Lamps window appears.
2. Make sure that the lamps you want to remove are switched off:  
The **On** buttons must *not* be green. If necessary, click on the respective **On** buttons.
3. Close the Align Lamps window.
4. Unplug and remove the lamps. For EDLs, remove the coding plugs as well.

### **Tips for installing lamps**

- Use only Perkin Elmer lamps. You can use both hollow cathode and electrodeless discharge lamps.
- Do not touch the front window of the lamp; perspiration or other contamination can reduce the intensity of the radiation.
- If you install more than one lamp containing a particular element, for example, a single-element copper lamp and a multi-element lamp containing copper, the system will use the lamp in the holder with the lower number when you perform determinations of that element. Make sure that you place the preferred lamp in the holder with the lower number.
- With EDLs, connect each lamp plug to the socket that has the same number as the holder of the lamp.
- With EDLs, make sure that you use the correct coding plug and connect it to the receptacle on the lamp holder containing the EDL.
- If you use the wrong coding plug or connect the coding plug to the wrong receptacle, you will cause the spectrometer to malfunction and may cause irreparable damage to the lamp.

## **Installing sample trays**

A sample tray can be inserted into one of three locations in front of the spectrometer – on the left-hand side, in front of the atomizer compartment, and on the right-hand side. You can leave a tray permanently installed on the left-hand side if you wish. You cannot install a tray in front of the atomizer compartment or on the right-hand side as long as the furnace autosampler is installed. When the autosampler is swivelled to the standby position at the right, you can install a tray in front of the atomizer compartment.

The sample trays are robust and can support an accessory, such as a diluter, or a flame autosampler, or a flow injection system.

### **To install the tray**

1. Hold the tray with both hands and position it in front of the location into which you wish to insert it.
2. Insert the tray support posts into the guides at the base of the instrument and slide in the tray until it hits the stop. Then gently lift the front of the tray and slide it finally into place. You can feel it lock into place.

### **To remove the tray**

1. Gently lift the tray at the front and then pull it back to slide it out of the guides.

## Shutting down the system

**Note:** In an emergency, switch off the spectrometer using the operational on/off switch on the front panel.

The flame is automatically extinguished and the gases shut down in a safe sequence.

See page 4-14 for the procedure to follow after an emergency shutdown.

You must shut down each atomizer correctly before you shut down the spectrometer.

### To shut down the flame system

1. With the flame still burning, aspirate the correct rinsing solutions to rinse the burner system (see 'Extinguishing the flame' in the 'Burner System User's Guide' for detailed information).  
Either aspirate the solutions manually or use the flame autosampler.
2. Extinguish the flame and bleed the gas lines:
  1. In the **Flame Control** window, click on the **Flame** icon.  
The system uses a predetermined shutdown sequence to extinguish the flame safely.
  2. Shut down the gases to the spectrometer at source.
  3. In the **Flame Control** window, click on **Bleed Gases** to depressurize the gas lines.

### To shut down the furnace system

1. Flush the furnace autosampler rinsing system (see 'Autosampler maintenance procedures' in the 'Graphite Furnace User's Guide' for detailed information).
2. Shut down the furnace gases to the spectrometer at source.

### To shut down the spectrometer

1. Exit AA WinLab:  
In the **File** menu, click on **Exit**.
2. Switch off the spectrometer at the operational on/off switch on the front panel.
3. Switch off any accessories.

4. If you do not want to use the computer for other tasks, shut down the computer and printer as described in their user's guides.
5. Empty the burner drain vessel (see 'Emptying the drain vessel' in the 'Burner System User's Guide').
6. Empty the furnace autosampler waste bottle.

**Note:** *Dispose of hazardous or corrosive solutions properly and refer to your local safety regulations for proper disposal procedures.*

## **Emergency shutdown**

If a situation arises where you have to shut down the system quickly and you do not have time to carry out the correct procedure, you can use the procedure described below.

*Note: This procedure is for use in an emergency only. Do not use it as a routine shutdown procedure.*

1. Switch off the spectrometer at the operational on/off switch on the front panel.
2. Shut down the gases to the spectrometer at source.

When you return to the spectrometer, proceed as follows to shut down the system properly.

3. If you were previously running concentrated salt solutions, organic solvents, or solutions containing cyanide, copper, silver, or mercury:
  1. Remove the burner head (*see the 'Burner System User's Guide'*).
  2. Pour a large quantity of deionized water slowly through the neck of the spray chamber to thoroughly flush the chamber and the drain system.
4. Check the burner head.  
If there are deposits along the slot, clean the burner head (*see the 'Burner System User's Guide'*).
5. Reinstall the burner head.
6. Switch on and set up the spectrometer system.
7. Ignite the flame.
8. Extinguish the flame properly.

# ***Spectrometer: System Maintenance***

---

# **5**

<b><i>Contents</i></b>	<b><i>page</i></b>
Perkin Elmer Service .....	5-3
Troubleshooting .....	5-3
Maintenance checklist .....	5-4
Spectrometer maintenance procedures .....	5-5
Cleaning the windows of the optical system .....	5-5
Maintaining the electrical components .....	5-7
Changing the air filters .....	5-9
Replacing the deuterium lamp .....	5-10

## Perkin Elmer Service

**Warning: Unauthorized Adjustments and Servicing**

W1.2

Only a Perkin Elmer service engineer or similarly trained and authorized person should be permitted to service the instrument.

- Do not attempt to make adjustments, replacements, repairs, or modifications to this instrument except as described in the documentation supplied with the instrument.

**If the spectrometer system does not function correctly...**

- Switch off the spectrometer system and make sure that it cannot be operated.
- Contact your local Perkin Elmer office.

You should only perform the maintenance procedures described in this guide. For any other maintenance or service contact your local Perkin Elmer office to arrange for a service engineer to visit.

In particular you should only allow a Perkin Elmer service engineer or similarly trained and authorized person to perform any work on:

- electrical components inside the instrument;
- the gas connections and gas system pneumatics;
- the monochromator and all other optical system components.

**Before the service engineer arrives:**

1. Make sure that the instrument and work area are clean.
2. Empty corrosive solutions from the burner drain vessel.  
Then add water to the drain trap to activate the function.
3. Empty the furnace autosampler waste bottle.

## Troubleshooting

See the *Troubleshooting Guide* supplied with the AA WinLab software. This guide includes information on problems that can arise with both the software and the spectrometer system.

## Maintenance checklist

<b>General preventative maintenance procedures</b>	
Instrument and work area	For safety reasons, and to avoid contaminating new samples, you should make sure that the instrument and the work area are always absolutely clean. This is especially important when you are working with toxic substances and when you are measuring trace amounts of any elements. Wipe up spills immediately before they can cause further contamination or damage. Before you leave the instrument standing idle for any period of time, make sure that all the components are clean.
Fume ventilation	Make sure that the main fume ventilation system functions correctly. Always switch it on when you are using the instrument.
Environmental conditions	Make sure that the environmental conditions are as described in ' <i>Laboratory Requirements</i> '.

<b>Spectrometer, maintenance checklist</b>	
External surfaces	Wipe over the surfaces with a lint-free cloth moistened with a dilute solution of laboratory detergent. The surfaces are resistant to dilute acids and alkalis, and to a lesser extent to strong acids and alkalis and organic solvents.
Air filters	Fit new filters yearly or more often in a dusty environment ( <i>procedure: page 5-9</i> ).
Atomizer compartment windows, lamp end windows	These windows do not normally require cleaning but if they become marked, e.g., with fingerprints, you can clean them ( <i>procedure: page 5-5</i> ).

## Spectrometer maintenance procedures

### **Procedures described**

- Cleaning the windows.
- Changing the air filter.
- Resetting the circuit breaker and changing the fuses.
- Replacing the deuterium lamp.

### **Cleaning the windows of the optical system**

---

#### **Caution: Risk of damage to optical components**

*Never touch or clean the surfaces of the grating or mirrors.*

*If mirror surfaces are scratched they must be resurfaced.*

*Do not allow cleaning fluids to come into contact with any of the reflecting surfaces.*

---

Most of the optical components are under the main cover of the spectrometer and are protected by the black optics covers.

Usually these components require no maintenance, provided the laboratory fume ventilation system is working correctly and the environmental conditions are as described in 'Laboratory Requirements'.

The only optical parts that are exposed are:

- The atomizer compartment windows.
- The windows of the lamps.
- The furnace windows. Cleaning the furnace windows is described under *System Maintenance* in the *Graphite Furnace User's Guide*.

These windows are important parts of the whole optical system. Handle them carefully, as you would any sensitive optical component. Do not directly touch the windows.

If mirror surfaces need cleaning, this must be done by a Perkin Elmer service engineer or a similarly authorized and trained person who knows the special techniques required.

**To clean the windows:**

This must be done only by a skilled person who knows the special techniques required.

*Note: The more you rub the surface of the windows, the more chance there is of scratching them. You must change scratched windows.*

1. Blow dust off optical surfaces with clean dry air before you attempt to clean them. Do not rub the surfaces.
2. Clean the window surfaces with a soft lint-free cloth moistened with spectroscopic grade alcohol. The window must dry quickly to avoid leaving drying marks.

**Keep the cleaning solutions and water away from other parts of the instrument.**

## Maintaining the electrical components

**Warning: Electrical Hazard**

W2.3

- To prevent potential injury to yourself and damage to the instrument, switch OFF all instruments in the system and disconnect them from the line power supply before you perform the procedures described below.

### Parts that may be serviced by the user

The only parts that you should attempt to maintain are those described below. The electronic components in the instrument do not contain any parts which you can maintain yourself. If there is a problem, contact your local Perkin Elmer office.

If any fuses blow frequently or if the automatic circuit breaker switch is activated frequently, contact your local Perkin Elmer office for advice.

### Resetting the automatic circuit breaker

The spectrometer has an automatic circuit breaker which prevents the system from drawing too much current.

When the circuit breaker is activated, the switch flips to the off position (marked O). This can happen while an analysis is running.

1. To reset the switch, wait a few seconds then put the switch to the on position (marked I).

If this switch is activated frequently, contact your local Perkin Elmer office for advice.

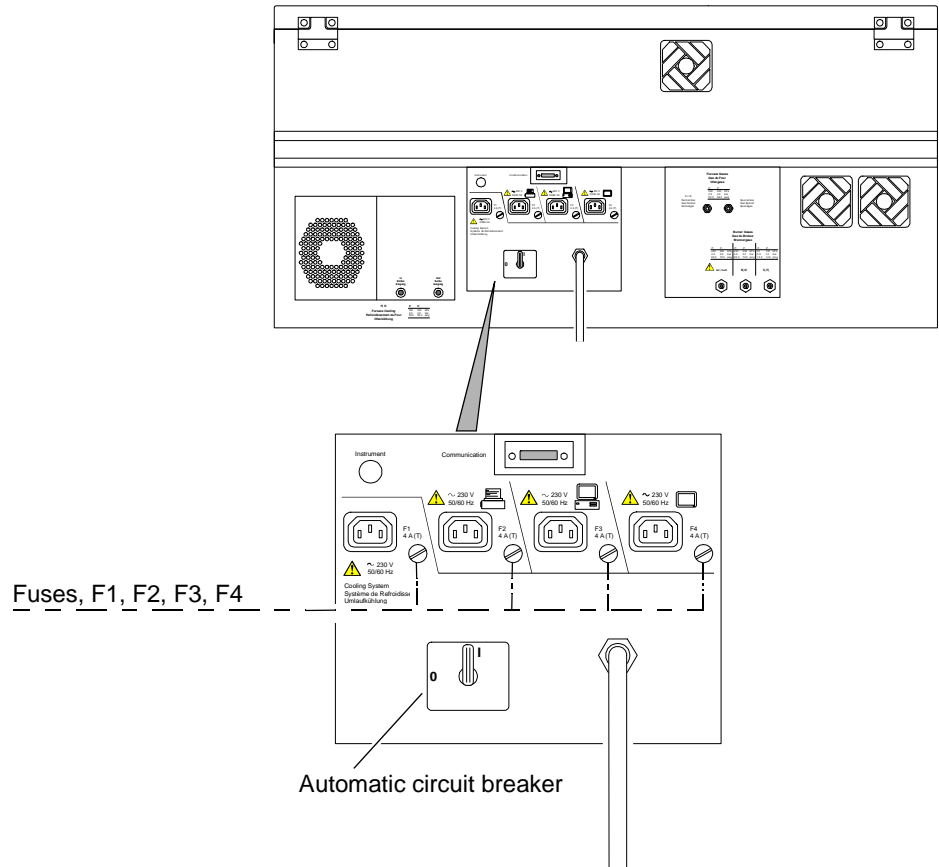
### Changing the fuses

**Warning: Electrical Hazard**

W2.9

- Use only fuses with the correct current rating and of the specified type for replacement.
- Do not use makeshift fuses and do not short-circuit the fuse holders.

1. Switch off the spectrometer and all other instruments in the system as described in the relevant guides.
2. Switch the circuit breaker on the rear panel of the spectrometer to the off position (marked O).

3. Disconnect the spectrometer's line power cord from the line power supply.
4. Disconnect all the line power cords from the sockets at the rear of the spectrometer.
5. To change the fuses numbered F1, F2, F3, F4:
  1. Insert the tip of a suitable screwdriver in the slit of the fuse holder and unscrew the fuse holder.
  2. Remove the blown fuse, then insert a new one of the correct type and rating. The correct rating is shown on the label of the fuse on the rear panel of the spectrometer.
  3. Screw the fuse holder into the correct socket.
6. Reconnect the line power cords correctly.


## Changing the air filters

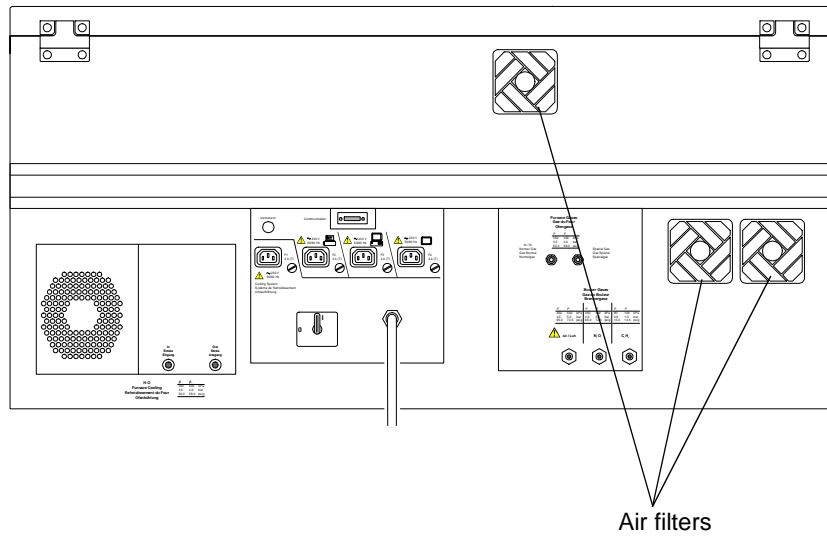To make sure that the electrical components in the spectrometer are properly cooled, the air filters must allow sufficient air into the instrument. Check the filters regularly and change them if they are dirty.

You should not need to change the filters more than once every year if the environmental conditions are as described under '*Laboratory Requirements*'. Replacement filters are available; see '*Parts and Supplies*' for the part number.

1. Carefully pull off the snap-on plastic grid that holds the air filter in place.
2. Replace the filter and push the grid back into place.

## Replacing the deuterium lamp

The deuterium lamp for the background corrector has a limited lifetime. After a prolonged period of operation the lamp will exhibit decreasing radiant intensity and will ultimately fail.

The lifetime of the deuterium lamp will be prolonged if you do not switch it on and off frequently.

You can replace the deuterium lamp assembly yourself. The replacement lamp is fitted in a prealigned mount.

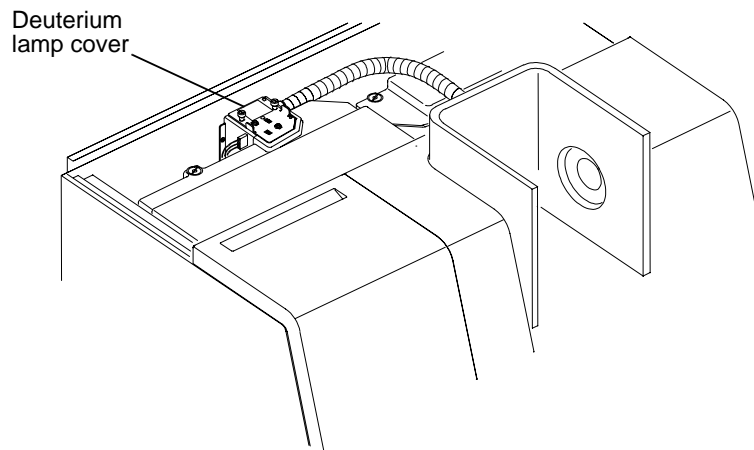
Figure 5-1. Location of Deuterium Lamp inside the Spectrometer


**Warning: Electrical Hazard**

W2.3

- To prevent potential injury to yourself and damage to the instrument, switch OFF all instruments in the system and disconnect them from the line power supply before you perform the procedures described below.

1. Switch off the spectrometer at the operational on/off switch on the front panel.
2. Switch the circuit breaker on the rear panel of the spectrometer to the off position (marked O).
3. Disconnect the spectrometer from the line power supply.

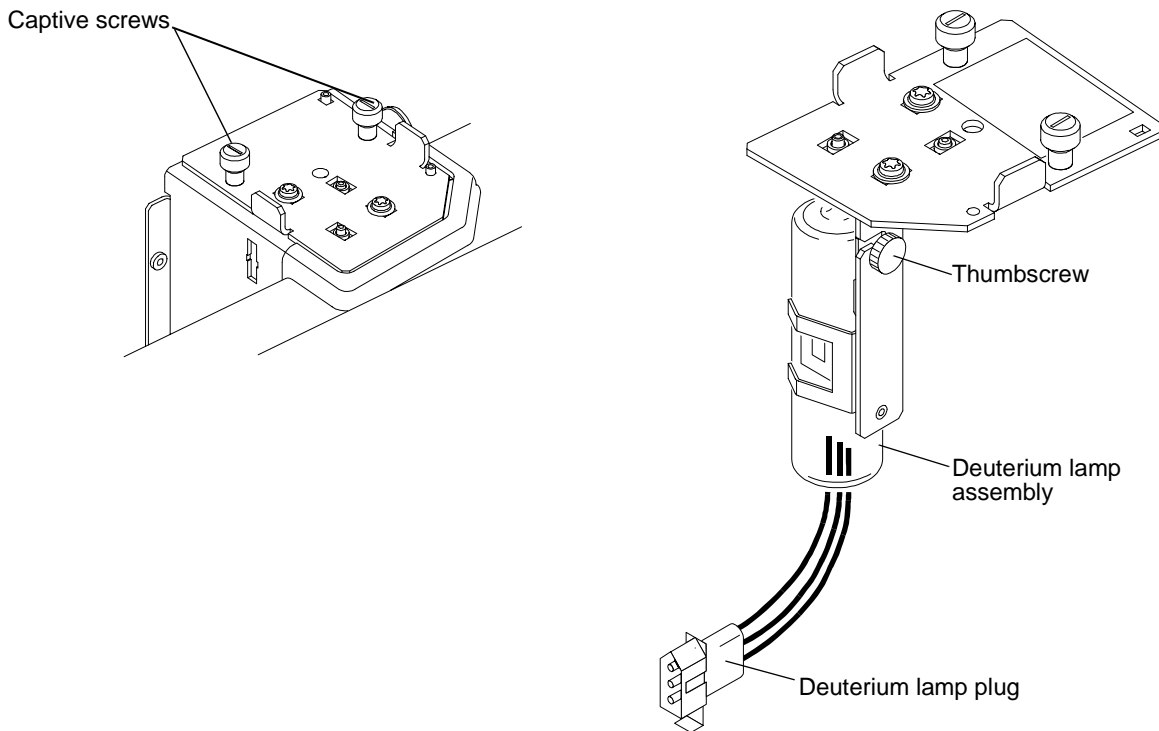
*Figure 5-2. Deuterium Lamp Cover and Bracket Assembly*

4. Open the top cover of the spectrometer:
  1. Remove the two securing screws; one on the left and one on the right.
  2. Swing back the cover to the open position.
5. Remove the deuterium lamp assembly:
  1. Undo the two captive screws on the deuterium lamp cover.
  2. Carefully lift up the deuterium lamp cover.  
The deuterium lamp assembly is attached to a bracket underneath the cover.
  3. Carefully place the complete assembly on top of the black optics cover.
  4. Pull out the deuterium lamp plug from the socket in the lamp compartment.  
**Do not pull on the deuterium lamp leads.**
6. Slacken the thumbscrew on the bracket and remove the old lamp assembly.  
Save the thumbscrew for use with the new lamp assembly.

7. Unpack the new lamp assembly, taking care to hold it only by the metal mount to prevent fingermarks on the lamp envelope.
8. Slip the slot at the base of the lamp mount over the stud on the lamp bracket and secure the lamp assembly with the thumbscrew saved in step 6.
9. Wipe the lamp envelope with a soft cloth moistened with alcohol to remove dirt, since this would otherwise be burned in when the lamp is hot.
10. Insert the deuterium lamp plug into the socket in the lamp compartment. The plug is asymmetric and can only be inserted in one direction.
11. Reinstall the lamp assembly:
  1. Carefully lower the deuterium lamp assembly into the lamp compartment until the lamp cover fits in place.
  2. Screw in the two captive screws on the lamp cover.
12. Close the top cover:
  1. Swing down the cover to the closed position.
  2. Screw in the two securing screws; one on the left and one on the right.

Since the deuterium lamp is fitted with a prealigned mount, you do not normally need to align the lamp.

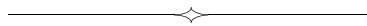**Note:** Do not attempt to make adjustments to the lamp bracket assembly. If there is indication that the alignment is not correct (e.g., low energy for the BG bar graph in the Align Lamps or Continuous Graphics windows), please contact your Perkin Elmer office.

# Parts and Supplies

---

# 6

<b>Contents</b>	<b>page</b>
Parts provided . . . . .	6-3
Replacement parts . . . . .	6-5
Obtaining supplies, replacement parts, and accessories . . . . .	6-5

**Parts provided**

Perkin Elmer reserves the right to alter the schedule of parts provided without prior notice.

<b>Qty.</b>	<b>Item</b>	<b>Part No.</b>
1	AAAnalyst 800 atomic absorption spectrometer. . . . .	B312-0081
1	Read cable. . . . .	B312-1302
3	Special line cords . . . . .	B310-0003
1 pack	Replacement air filters . . . . .	B050-2706
4	Plugs, for handle screw holes . . . . .	B015-3803
1 pack	Replacement fuses . . . . .	B312-1303
1	AS-800 furnace autosampler . . . . .	B300-1500
1	Swing-arm mount for autosampler. . . . .	B312-0923
	<i>Further parts provided with the furnace autosampler are listed under 'Parts and Supplies' in the 'Graphite Furnace User's Guide'.</i>	
1	Cooling system . . . . .	B019-8072
1	Fuel hose assembly (red) . . . . .	B015-4984
1	Air hose assembly (black) . . . . .	B015-4987
1	Nitrous oxide hose assembly (blue) . . . . .	B015-4990
10 m	Special gas tube (FLEXO) . . . . .	B050-8264
1	Y-connector. . . . .	0992-0179
1	Drain vessel . . . . .	B312-1306
1	Sample tube . . . . .	B015-6272
5	Corkpreen gaskets for burner end cap . . . . .	0047-2014
1	O-ring for burner end cap . . . . .	B015-8690
1	O-ring for the spray chamber . . . . .	B004-7207
1	Burner head cleaning tool. . . . .	0303-1573
1 pack	Graphite tubes. . . . .	B300-0409
1 set	Graphite contacts . . . . .	B300-0410

<b>Qty.</b>	<b>Item</b>	<b>Part No.</b>
1 pack	Cleaning swabs . . . . .	B050-5368
1	Grips for graphite tube . . . . .	0990-7925
1 set	Tools, for changing contacts and furnace windows . . .	B019-3347
1	Instrument mirror . . . . .	B085-1900
1	Replacement O-ring, for front contact housing and optical sensor in the rear contact housing . . . . .	B013-2870
1	Documentation package	B312-1326

## Replacement parts

### Obtaining supplies, replacement parts, and accessories

Supplies, replacement parts, and accessories can be ordered directly from Perkin Elmer, using the part numbers quoted in this guide.

Perkin Elmer's catalog service offers a full selection of high-quality atomic spectroscopy supplies through the *Atomic Spectroscopy Supplies Catalog*.

To place an order for supplies and many replacement parts, request a free catalog, or ask for information:

- If you are located *within the U.S.*, call toll free **1-800-762-4002**, 8 a.m. to 8 p.m. EST. Your order will be shipped promptly, usually within 24 hours.
- If you are located *outside of the U.S.*, call your local Perkin Elmer sales or service office.

### What is listed in this section

This section lists the most important replacement parts for your spectrometer system. Parts listed under *Parts provided* are not repeated in this section.

Information about other parts and accessories for the system is given in the current Perkin Elmer brochures and price lists.

### Background corrector

Description	Part No.
Deuterium lamp . . . . .	B050-7404

<b>Contents</b>	<b>page</b>
Introduction .....	7-3
Space requirements for the spectrometer system .....	7-4
Magnetic field .....	7-7
Electrical requirements .....	7-8
Environmental conditions .....	7-9
Fume ventilation .....	7-10
Gas supply requirements .....	7-11
Burner gases .....	7-12
Furnace gases .....	7-14
Cooling the spectrometer system .....	7-15
Lamps .....	7-16
Computer and printer .....	7-16

## **Introduction**

The Analyst 800 spectrometer has some special requirements that you must consider when setting up the laboratory. The laboratory must fulfil these requirements, described in this chapter, before you can operate the spectrometer correctly.

Brochure no. L-301 – *Preparing Your Laboratory for Perkin Elmer Atomic Spectroscopy Instruments* – also contains useful information. This brochure is available from Perkin Elmer.

## **Space requirements for the spectrometer system**

The illustrations on the following pages show the dimensions of the spectrometer and the space that you require to accommodate the spectrometer, computer, and ancillary equipment.

- Place the spectrometer on a sturdy bench or trolley capable of sustaining the weight. A movable table is very convenient, but make sure that the wheels can be securely locked.
- You should have sufficient room around the instruments to reach all the connections on the rear and reach the electricity and gas supply points.
- The laboratory ventilation system needs space above the spectrometer.

### **Computer and printer**

Place the computer near the spectrometer in a convenient position for working with both instruments. A movable table, specially designed for a computer and printer, may be most convenient.

The printer needs to be in a convenient place with enough room for the paper supply to stand and feed smoothly into the printer.

### **Accessories**

Make sure that there is sufficient room for any accessories near to the spectrometer.

For example, we recommend that you place the (optional) FIAS flow injection system on a movable table in front of the atomizer compartment to minimize the length of the sample transfer tube. Make sure that the wheels of the table can be securely locked.

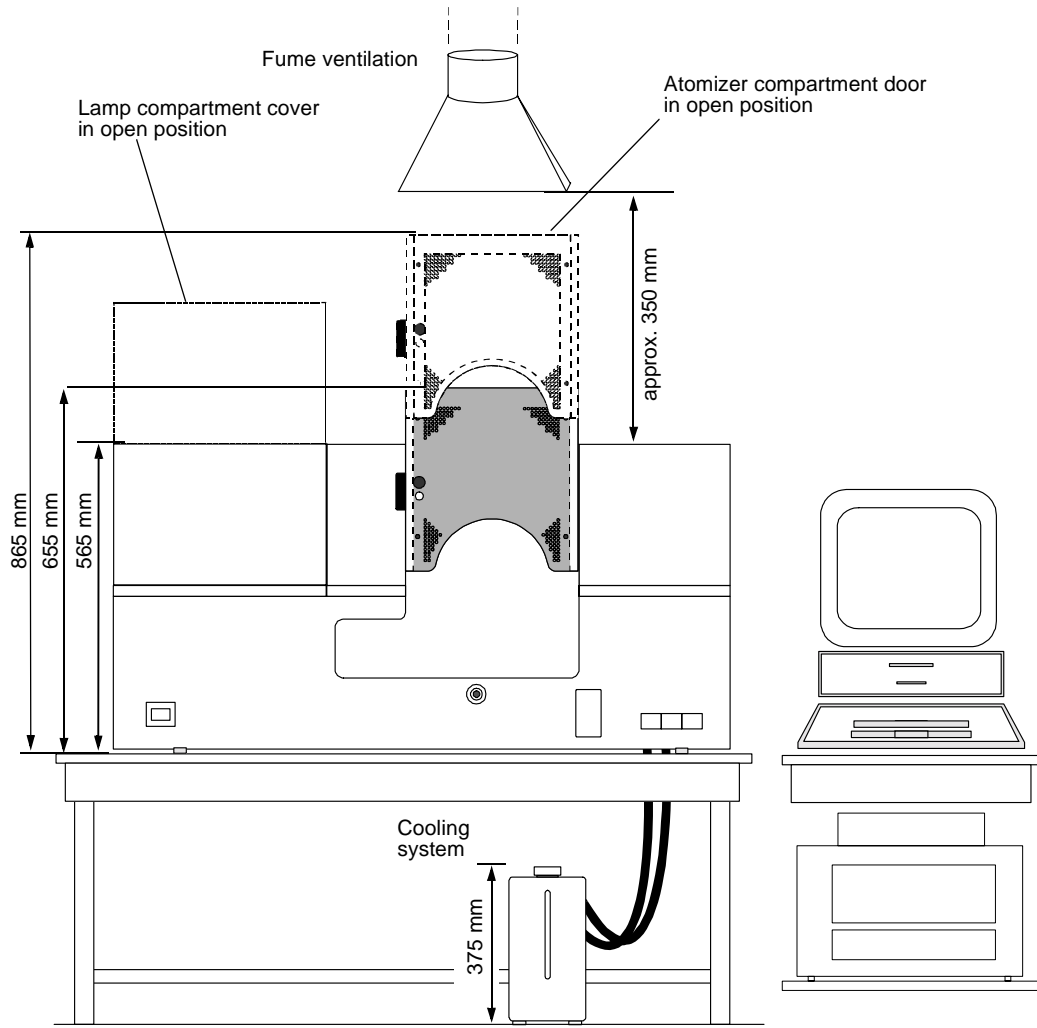
Figure 7-1. Space Requirements for the Spectrometer System (1)

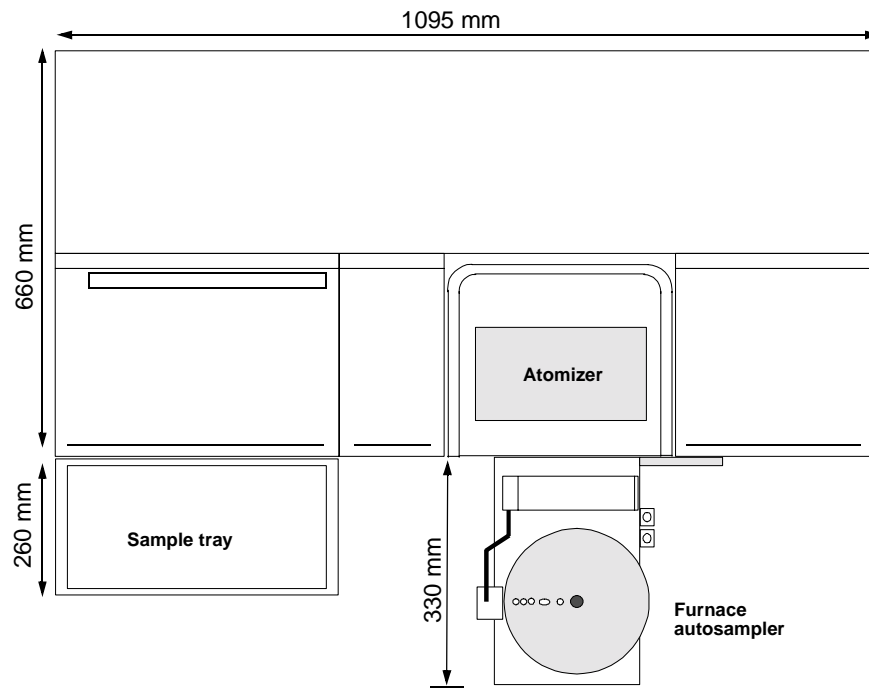
Figure 7-2. Space Requirements for the Spectrometer System (2)



## Magnetic field
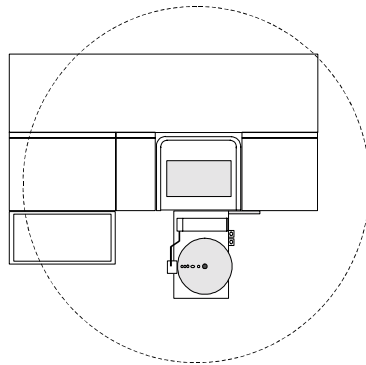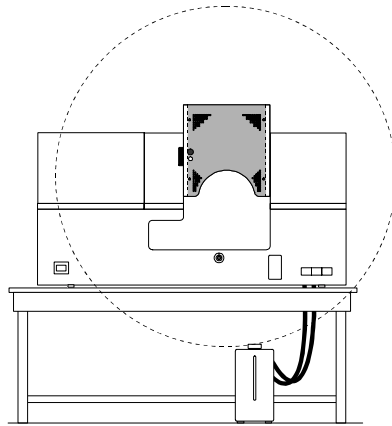
### Zeeman graphite furnace systems

The electromagnet generates a strong magnetic field inside the THGA furnace during measurement cycles.

There is a weak, accessible, stray magnetic field around the graphite furnace. This stray magnetic field can affect heart pacemakers and metallic implants up to 0.6 meter from the furnace in all directions. This stray field can also affect other instruments, watches, and magnetic storage devices.

- Do not operate the furnace when persons wearing electronic heart pacemakers, or having other metallic implants, are present.
- Anyone wearing a heart pacemaker or having metallic implants should remain at least 0.6 meter away from the furnace – in any direction – while the furnace is operating.

#### Strong Magnetic Field

Safety distance of 0.6 meter (24 inches) radius around the furnace magnet for persons wearing heart pacemakers

## **Electrical requirements**

Two versions of the AAnalyst 800 AA spectrometer are available to suit different line power supplies.

### **USA version – 240 V, 60 Hz**

The line power cable has a L6-20P, NEMA-configuration twistlock connector (3 wire – line, neutral, grounding), rated for 240 V/20 A, the cable is 3 meters long.

### **General version – 230 V, 50 Hz**

The line power cable has a 3-pole CENELEC connector (line + neutral + protective earth) rated for 32 A. The cable is 3 meters long.

### **General requirements**

The spectrometer requires a stable, 50 Hz or 60 Hz line power supply at the correct voltage and current rating that has a correctly wired protective earthing system (ground connection).

The line power supply should be free of line transients in excess of 50 V peak. If the electrical supply voltage produces large AC line voltage fluctuations, a qualified electrician should install a voltage regulator between the electrical outlet and the instrument.

Power consumption during furnace operation is approximately 5000 VA, at an isothermal furnace temperature of 2600 °C.

The line power supply point must be within 3 meters of the rear of the spectrometer.

We recommend that you connect the spectrometer, computer, printer, and any accessories to the same phase of the line power supply and the same protective earth.

The line power supply must conform with local safety regulations and be checked by a qualified electrician before you connect the instrument to line power.

See the guides for the computer and printer and any other instruments for information about their power requirements.

## Environmental conditions

**Warning: Explosive Atmosphere**

W1.3

*This instrument is not designed for operation in an explosive atmosphere.*

**General**

The location you choose for the spectrometer must fulfil the criteria listed below:

- Indoors.
- Free of dust, smoke and corrosive fumes.
- The work surface must be level, flat, clean, dry, and free from vibration. It should be strong and stiff enough to bear the weight of the spectrometer system over a long period without warping or bending.
- Out of direct sunlight.
- Away from radiators and heaters.
- Near to the electricity and gas supply points.
- Near to a source of cooling water for the graphite furnace.
- Have a suitable fume ventilation system.
- Have sufficient room to work comfortably with the instruments, and allow you to reach the connections at the rear of the instruments.

**Temperature**

- Between +15 °C and +35 °C.

**Humidity**

- Between 20% and 80% relative humidity (non-condensing).

## **Fume ventilation**

You must provide an efficient fume ventilation system to remove the fumes generated by the atomizer (flame, graphite furnace, or the QTA-cell of FIAS systems).

The specification for a suitable system is described in brochure no.

L-301 – *Preparing Your Laboratory for Perkin Elmer Atomic Spectroscopy Instruments.*

This brochure is available from Perkin Elmer.

**Warning: Toxic Gases – Fume Ventilation System**

*Without adequate ventilation, potentially toxic vapors can build up in the laboratory.*

- *Your laboratory must have a reliable fume ventilation system before you use this instrument.*

## Gas supply requirements

The permanent installation of gas supplies is the responsibility of the user and should conform to local safety and building regulations.

Perkin Elmer recommends that cylinders be stored outside the laboratory and that gases be piped to the spectrometer system in approved gas lines.

- The user must provide the gas supplies, regulators, connectors, and valves. A double-stage or heated regulator is recommended for nitrous oxide. Either the gas supplier or Perkin Elmer can supply the correct regulators and valves.
- You must be able to reach the on/off valves easily and see the pressure indicators.
- Consult the chapter entitled '*Safety Information*' for recommended safety procedures.

**Warning: Compressed Gases**

*High pressure gas cylinders can be dangerous if mishandled or misused.*

- *Always handle gas cylinders with caution and observe you local regulations for the safe handling of gas cylinders.*

## Burner gases

Two types of flame are commonly used for atomic absorption analyses with flame atomization: the air/acetylene flame and the nitrous oxide/acetylene flame. The specifications and requirements for the burner gases are given below:

Table 7-1. Specifications and Requirements of the Burner Gases

Gas	Purity	Outlet Gauge Pressure					
		kPa		bar		psig	
		$P_{\min}$	$P_{\max}$	$P_{\min}$	$P_{\max}$	$P_{\min}$	$P_{\max}$
Acetylene (ethyne, $C_2H_2$ )	99.996%, atomic absorption grade in acetone	90	100	0.9	1.0	13.0	14.5
Air	oxygen $\leq 20\%$ ; particle free; water free	450	500	4.5	5.0	65.0	72.5
Nitrous oxide (dinitrogen monoxide, $N_2O$ )	99.996%	450	500	4.5	5.0	65.0	72.5

- Set the outlet gauge pressure for each gas to a value just below the maximum value ( $P_{\max}$ ).
- Always set the same outlet gauge pressure for air and nitrous oxide.
- **Never** set the outlet pressure to a value exceeding the maximum pressure ( $P_{\max}$ ).

**Warning: Acetylene – Explosion Hazard**

W5.7

Acetylene can decompose explosively at pressures higher than 103 kPa/  
1.03 bar/15 psig.

- Always make sure that the acetylene outlet gauge pressure is below this value.

### **Air supply**

An air compressor is recommended for the air supply to the burner system. Cylinders of compressed air should only be used as an emergency or short-term solution. Since compressors are generally noisy, we recommend that you locate the compressor outside the laboratory.

#### **Warning: Flashback Hazard**

W5.9

*Perkin Elmer burner systems are designed for use with compressed air. The use of oxygen can cause an explosion in the burner system, and oxygen-enriched air can cause a flashback of the flame.*

- *Never use oxygen or oxygen-enriched air with Perkin Elmer burner systems. Use only compressed air.*

## Furnace gases

### Inert gas

The graphite furnace requires a supply of inert gas to prevent the tube and the analyte atoms from being oxidized when the tube is heated.

Argon is the recommended inert gas.

If you cannot obtain argon, you can use purified nitrogen, but nitrogen does not provide such efficient protection as argon at temperatures above 2000 °C.

Nitrogen can also react with the samples and the graphite tube at typical atomization temperatures.

#### Warning: Toxic Gases

W6.1

If you use nitrogen as the inert gas, toxic cyanogen gas (CN)<sub>2</sub> will be produced when the furnace temperature is above 2300 °C.

- Make sure that the main laboratory fume ventilation system removes all gases and fumes emitted by the furnace.

### Special gas

For some applications you may want to use a special gas. A typical example is the use of air during one of the pyrolysis steps to better ash organic materials.

Table 7-2. Specifications and Requirements of the Furnace Gases

Gas	Purity	Outlet Gauge Pressure						Max. Flowrate
		kPa		bar		psig		L/min
		$P_{\min}$	$P_{\max}$	$P_{\min}$	$P_{\max}$	$P_{\min}$	$P_{\max}$	1.2
Argon	99.996%	350	400	3.5	4.0	50.0	58.0	1.2
Nitrogen	99.996%	350	400	3.5	4.0	50.0	58.0	1.2
Special	particle free; water free	350	400	3.5	4.0	50.0	58.0	1.2

- Set the outlet gauge pressure for each gas to a value just below the maximum value ( $P_{\max}$ ).
- **Never** set the outlet pressure to a value exceeding the maximum pressure ( $P_{\max}$ ).

## **Cooling the spectrometer system**

### ***Furnace cooling***

The graphite furnace uses water cooling to allow the graphite tube to return to its equilibrium temperature (approximately ambient) quickly between measurements (*see 'Cooling the furnace' in the 'Graphite Furnace User's Guide'*).

### ***Spectrometer cooling***

Cooling coils are incorporated in the spectrometer's power supply to maintain the electronic components at a comfortable working temperature.

### ***Cooling system***

Perkin Elmer offers a recirculating cooling system; it is included with certain spectrometer systems and is offered as an option for others. It provides coolant at a constant temperature and flow, to allow the most reproducible analytical conditions. It also makes the system independent of external water supplies and saves water (*see 'Cooling the furnace' in the 'Graphite Furnace User's Guide'*).

Power requirements: 230 V; 50/60 Hz (electrical power is supplied by the spectrometer)

Part Number: B019-8072

## **Lamps**

These are often called primary source lamps, atomic absorption source lamps, spectroscopic lamps, or line sources.

You need to provide the lamps for your particular applications.

Brochure no. L-656 – *Hollow Cathode Lamps and Electrodeless Discharge Lamps for Atomic Absorption*, contains descriptions of all the lamps that are available.

## **Computer and printer**

Perkin Elmer offers a suitable computer for controlling the spectrometer and a suitable printer. Contact your local Perkin Elmer office for more information. If you are not using the computer offered with the spectrometer by Perkin Elmer, or an identical model with the correct specification, Perkin Elmer cannot accept responsibility for incorrect functioning of the spectrometer system.

# ***Spectrometer: System Description***

---

# **8**

<b><i>Contents</i></b>	<b><i>page</i></b>
The AAnalyst 800 Spectrometer System .....	8-3
The optical system .....	8-6
The gas and cooling connectors .....	8-8
The electrical connectors and switches .....	8-10
Technical data .....	8-13

## **The AAnalyst 800 Spectrometer System**

### **Concept**

The AAnalyst 800 spectrometer from Perkin Elmer is a compact, high performance atomic absorption spectrometer incorporating a burner system for flame atomization and a graphite furnace for electrothermal atomization. The spectrometer is capable of fully automatic single-element analyses. The complete system is operated from the associated computer using the AA WinLab software.

### **Features of the AAnalyst 800 System**

- Two fully integrated atomizers: a burner system for flame atomization and a graphite furnace for electrothermal atomization. Atomizer changeover is automatic at the click of a mouse button.
- Transversely-heated graphite furnace featuring full Stabilized Temperature Platform Furnace (STPF) technology.
- Longitudinal Zeeman-effect background correction for electrothermal atomization.
- Deuterium background correction for flame atomization.
- Automated self-optimization of the burner system.
- High capacity furnace autosampler.
- Easy-to-use AA WinLab software with which you control the spectrometer system. AA WinLab also incorporates a range of powerful quality control functions.

### **System components**

The AAnalyst 800 is a compact benchtop spectrometer. The major components of the system are:

- The spectrometer, containing the optical system, the electronics, and the atomizer compartment.
- The burner system, mounted on a carriage that can be driven into and out of the atomizer compartment by software command (*see 'Burner system design' in the 'Burner System User's Guide' for more information*).
- The graphite furnace, mounted on a carriage that can be driven into and out of the atomizer compartment by software command (*see 'Furnace design' in the 'Graphite Furnace User's Guide' for more information*).
- The furnace autosampler, mounted on a swing arm in front of the atomizer compartment (*see 'Autosampler design' in the 'Graphite Furnace User's Guide' for more information*).
- The furnace cooling system (*see 'Cooling the furnace' in the 'Graphite Furnace User's Guide' for more information*).
- The computer and the associated AA WinLab software.
- The printer; this is optional but strongly recommended.

### ***System control***

To control the spectrometer and perform your analyses, you use the AA WinLab application program, which runs under the Microsoft Windows 95 operating system. The computer keyboard, the screen, and the mouse form the user interface.

When you switch on the spectrometer system it performs an initialization procedure to check the instrument. While the spectrometer is operating, the control system continually monitors many instrument functions. It generates appropriate messages when a problem occurs and prevents potentially hazardous situations from arising.

You can obtain more information about the application program in the AA WinLab online help.

### ***Accessories***

Perkin Elmer offers autosamplers for operation with the burner system, flow injection systems, diluters, and much more. Please contact your local Perkin Elmer office to obtain more information.

## **The optical system**

The AAAnalyst 800 uses a double-beam optical system. Front-surfaced, reflecting optical components are used throughout. The mirror surfaces are coated with silica for durability.

The lamp compartment has positions for up to eight spectral sources. The radiation from the requested source lamp is reflected by a movable mirror into the optical path.

For flame atomization a beam splitter splits the radiation into a sample beam and a reference beam. The sample beam passes through the atomizer while the reference beam passes around the atomizer and is conducted by a fiber optic. The sample and reference beams are measured simultaneously.

For electrothermal atomization the beam splitter is moved out of the optical path and a single-path optical configuration is used. Due to the Zeeman principle, the double-beam function is retained.

The Littrow-configuration monochromator has a large area reflecting grating blazed at two wavelengths; 236 nm for operation in the UV spectral range and 597 nm for operation in the visible spectral range. The monochromator offers a choice of three spectral slit widths. Two slit heights are available for each slit width. The high slits are used with the flame technique while the low slits are used with the graphite furnace technique. The detector is a segmented array of discrete photodiodes.

For continuum source background correction with the flame technique, the radiation from the deuterium lamp is merged with the source radiation at the beam splitter.

The inverse longitudinal Zeeman effect is used for background correction with the graphite furnace technique. This system of background correction is described in the '*Graphite Furnace User's Guide*'.

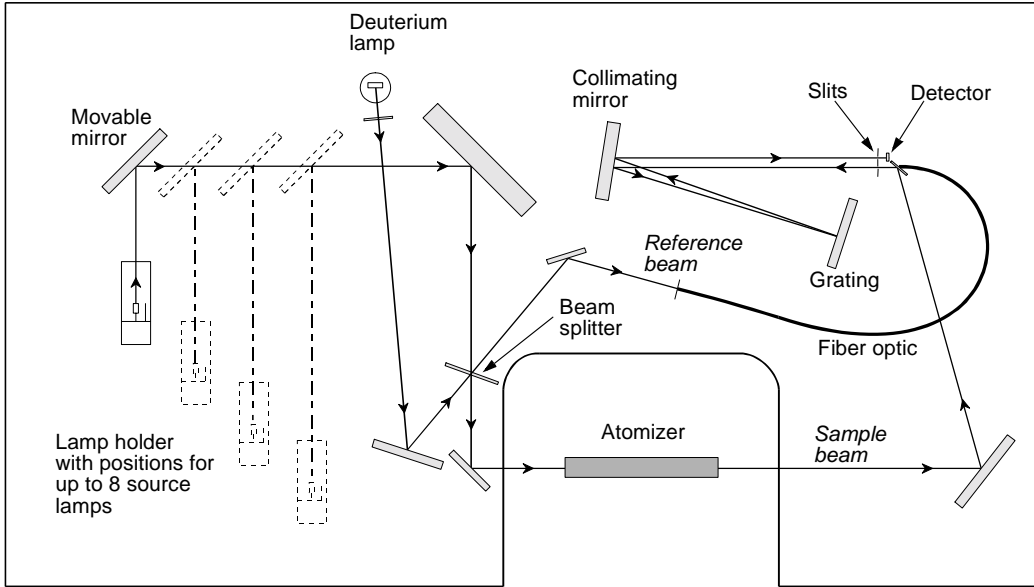
Figure 8-1. Optical System

**The gas and cooling connectors**

<b>Item</b>	<b>Designation</b>	<b>Function</b>
<b>Furnace Cooling</b>		
1	In	Fitting for connection of the cooling water inlet hose to the spectrometer.
2	Out	Fitting for connection of the cooling water outlet hose from the spectrometer.
<b>Furnace Gases</b>		
3	Ar/N <sub>2</sub> Normal Gas	Connector for the connection of the normal gas supply to the furnace (argon or nitrogen).
4	Special Gas	Connector for the connection of the special gas supply to the furnace. If you are not using a special gas, you must also connect the normal gas to this connector.
<b>Burner Gases</b>		
5	Air/Luft	Fitting for connection of the air supply to the spectrometer. The air supply is required for the burner and also for the pneumatic system in the spectrometer.
6	N <sub>2</sub> O	Fitting for connection of the nitrous oxide supply to the burner.
7	C <sub>2</sub> H <sub>2</sub>	Fitting for connection of the acetylene supply to the burner.

---

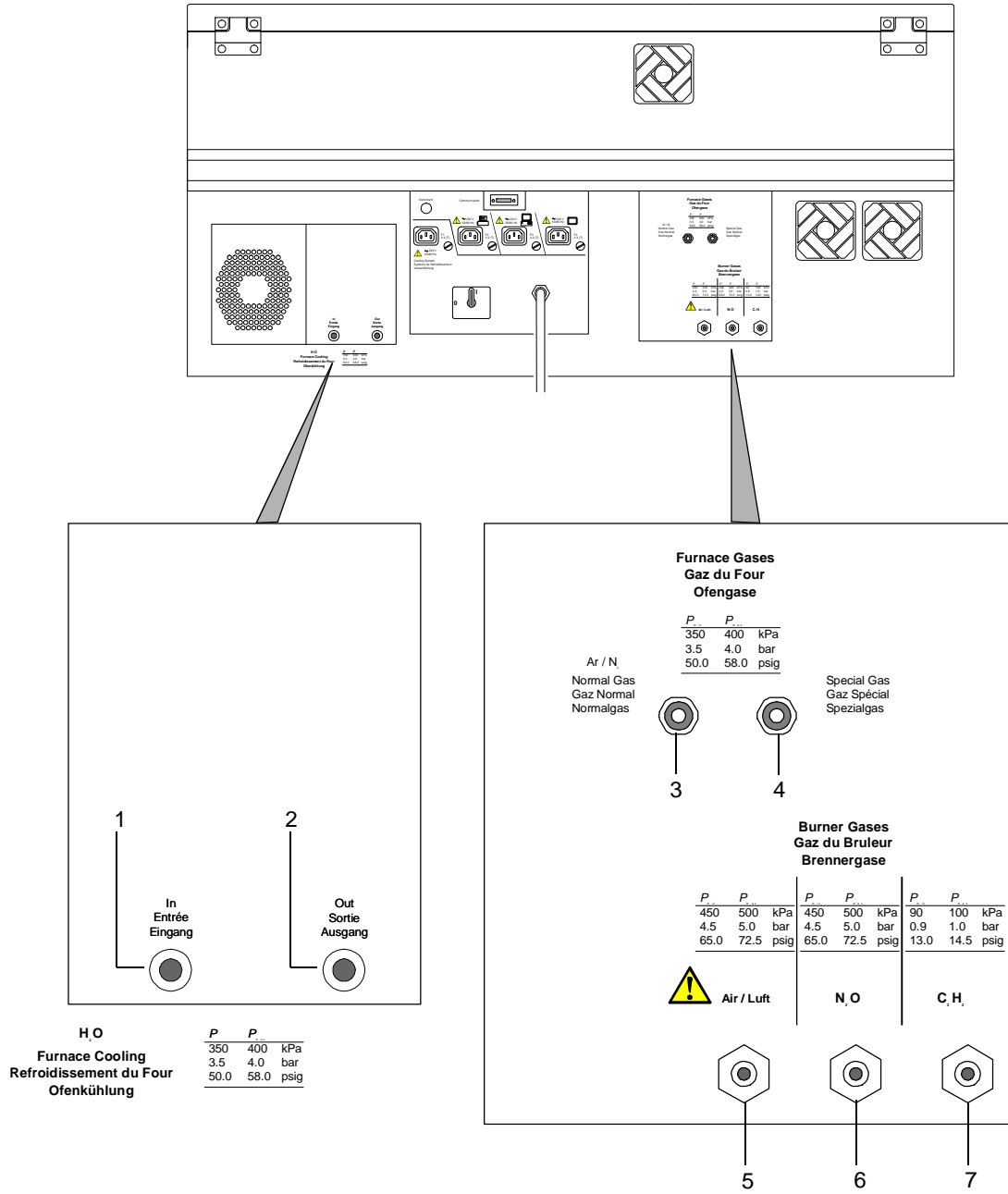
Figure 8-2. Gas and Cooling Connectors on the Rear Panel

## The electrical connectors and switches

### Connectors and controls on the rear panel

Item	Designation	Function
1	<b>Communication</b>	Socket for the communication cable connecting the spectrometer to the computer.
2	<b>Instrument</b>	Socket for a read cable to an external instrument to activate a read cycle.
3	<b>Cooling System</b> ~230 V 50/60 Hz	Power outlet for the cooling system. Connect the cooling system to this socket. <b>Do not connect other equipment to this power outlet.</b>
4	 ~230 V 50/60 Hz	Power outlet for a printer.
5	 ~230 V 50/60 Hz	Power outlet for a computer.
6	 ~230 V 50/60 Hz	Power outlet for a monitor.
7	○ 	Circuit breaker. An automatic switch that prevents the system from drawing too much current. It can also be used to isolate the spectrometer system from the line power supply.
8	---	Captive line power cord for the complete system.
9	<b>F1</b>	Fuse for the cooling system power outlet; 4 A slow-blow (T).
10	<b>F2</b>	Fuse for the printer power outlet; 4 A slow-blow (T).
11	<b>F3</b>	Fuse for the computer power outlet; 4 A slow-blow (T).
12	<b>F4</b>	Fuse for the monitor power outlet; 4 A slow-blow (T).

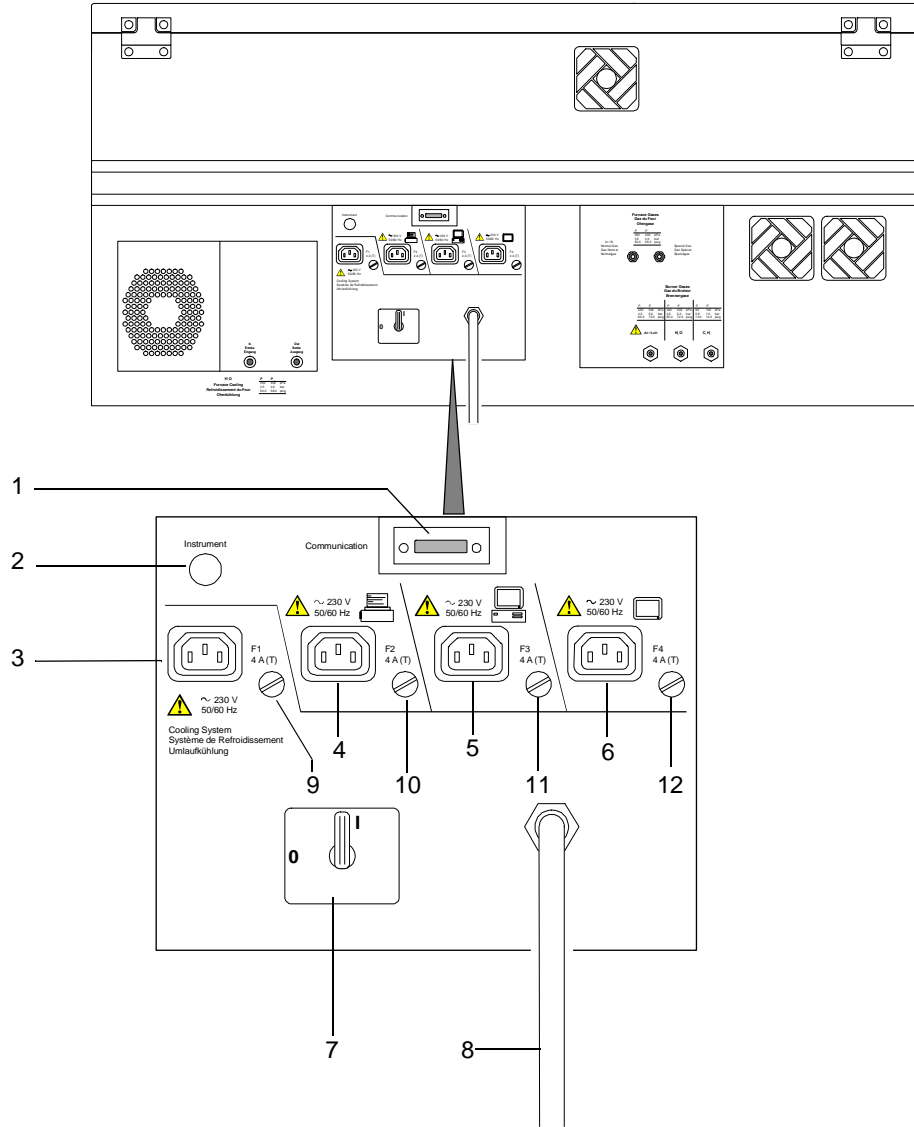
Figure 8-3. Electrical Connectors and Switches on the Rear Panel

### Switches and ports on the front panel

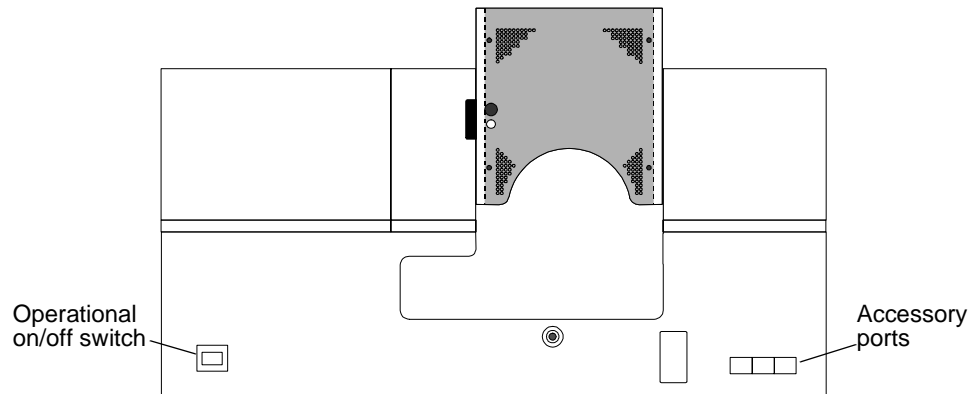
Figure 8-4. Switches and Ports on the Front Panel

### Operational on/off switches

#### Spectrometer

You normally switch the system on and off with the operational on/off switch on the front panel of the spectrometer.

#### Cooling system

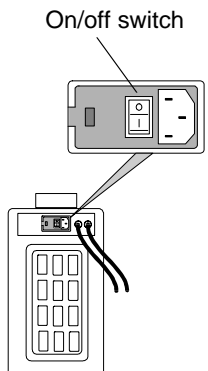
The cooling system requires a 230 V line power supply, which is supplied and controlled from the spectrometer. A special line power cord is provided for this purpose. Do not connect the cooling system to an external power source. You can leave the on/off switch on the cooling system in the 'on' position since the system is controlled from the spectrometer.

#### Accessory ports

Three accessory ports are located on the front panel of the spectrometer behind slide-up covers. You use these ports to connect certain accessories to the spectrometer, such as the furnace autosampler or a diluter.

The ports provide the electrical supply and control signals for operation of the accessories.

#### Cooling System

**Technical data**

<b>Principle</b>	Atomic absorption spectrometer. Computer controlled, using a special application program running under a graphical user interface. Real-time double-beam, all-reflecting optical system (single-path for furnace operation). Front-surfaced optics protected by a silica coating. AC modulated spectral line sources.
<b>Monochromator</b>	Littrow configuration grating monochromator. Grating: ruled area: 64 mm x 72 mm, 1800 lines/mm blaze wavelengths: 236 nm and 597 nm Focal length: 267 mm Slit widths 0.2, 0.7 and 2.0 mm Wavelength range: 190–870 nm Reciprocal linear dispersion: 1.6 nm/mm (nominal)
<b>Detector</b>	Segmented discrete solid-state detector with CMOS charge amplifier array; software-controlled automatic gain control (AGC).
<b>Lamps</b>	Line radiation sources. Positions for up to 8 lamps in lamp compartment. Single- and multi-element hollow cathode lamps (HCLs) and up to 4 electrodeless discharge lamps (EDLs) can be used.
<b>Background correction</b>	Continuum source (deuterium lamp) for the flame technique. Inverse longitudinal Zeeman effect for the graphite furnace technique.
<b>Atomizers</b>	Flame atomization: Premix burner system with pneumatic nebulizer and interchangeable burner heads; automatic changeover between air/acetylene and nitrous oxide/acetylene flames. Electrothermal atomization: Transversely-heated graphite furnace incorporating an electromagnet for longitudinal Zeeman-effect background correction.

<b>Signal handling</b>	<p>Time average, peak average, peak height, peak area.</p> <p>Measurement ranges:</p> <ul style="list-style-type: none"><li>absorbance: -0.5 to 2.000</li><li>concentration: 0.00001 to 9999 concentration units</li><li>expansion factor: 0.01 to 100</li><li>read time: 1 to 60 s</li><li>read delay: 0 to 60 s</li></ul>
<b>Calibration</b>	<p>Automatic zeroing, resloping, and recalibration.</p> <p>Calibration/recalibration using up to 15 calibration solutions ('standards'). Calibration coefficients determined by a least squares procedure.</p> <p>Analytical curve technique:</p> <ul style="list-style-type: none"><li>• Linear with zero or non-zero intercept.</li><li>• Non-linear, 2-coefficient, or 3-coefficient, zero or non-zero intercept.</li></ul> <p>Analyte addition technique:</p> <ul style="list-style-type: none"><li>• Zero or non-zero intercept.</li><li>• Additions calibrate (zero intercept).</li></ul>
<b>Computer</b> (minimum requirement)	<p>PC with Intel Pentium processor, 16 MB RAM, CD-ROM drive.</p> <p>Hard disk space: 30 MB minimum free for AA WinLab, Database and GPIB software, additional space required for storing results and methods.</p> <p>Microsoft Windows 95 operating system.</p> <p>Perkin Elmer offers a suitable PC for controlling the spectrometer system.</p>
<b>Application program</b>	<p>Special application program: AA WinLab.</p> <p>Online help function.</p> <p>Hard disk or floppy storage of measurement data, signals, and completed analytical methods.</p> <p>Special Reporter, Reformatter, and Library Manager applications for generating data reports and reformatting data for export to other data handling programs.</p>
<b>Graphics</b>	<p>High resolution display of the signals.</p> <p>Simultaneous display of the corrected analyte signal and the background signal.</p> <p>Calibration curve display.</p> <p>Storage and recall of transient absorbance signals.</p> <p>Hard copy printout of screen graphics on an associated printer.</p>
<b>Power requirements</b>	<p>230 V single-phase, 50 Hz, power consumption 5000 VA, -or- 240 V single-phase, 60 Hz, power consumption 5000 VA</p>

<b>Electrical protection</b>	As defined in EN 61010-1-1993 (IEC 1010-1): Insulation: Class I Insulation category (overvoltage category): II Pollution degree: 2
<b>Safety standards</b>	As defined in EN 61010-1-1993 (IEC 1010-1: 1990 + A1 + A2) and EN 61010-2-061:1996.  Compliant to the requirements of Canadian Standards Association (CSA) standards CAN/CSA-C22.2 No. 1010.1-92, Amendment 1 (1992), 2 (1995), and CAN/CSA 61010-2-061 (1995); File No. LR 45111.  Compliant to UL 3101-1 1993 and UL 3111-1 1994.  Certificate of Approval and GS mark from the LGA Bavaria, FRG, Institute of Equipment Safety Testing and Certification.
<b>EMC standards</b>	As defined in EN 61326-1:1997.
<b>Environmental requirements</b>	Ambient temperature: +15 °C to +35 °C, 20–80% relative humidity; non-condensing
<b>Dimensions</b>	Width: 1095 mm                      with handles: 1220 mm Height: 655 mm                      with door open: 865 mm Depth: 700 mm                      with handles: 770 mm with autosampler: 1040 mm
<b>Mass (weight)</b>	187 kg Spectrometer alone, not including the computer, autosampler, and cooling system.

# Translations

---

## Contents

	<i>page</i>
Symbols used on the instrument .....	T-3
Warning messages .....	T-11

## **Symbols used on the instrument**

Any of the following symbols can be used on the instrument:

<i>dansk</i>	<b><i>De markeringer, som benyttes på dette apparat</i></b> <i>Enhver af de følgende markeringer kan benyttes på apparatet:</i>
<i>Deutsch</i>	<b><i>Symbole am Gerät</i></b> <i>Eines der folgenden Symbole kann sich am Gerät befinden:</i>
<i>español</i>	<b><i>Pictogramas usados en el equipo</i></b> <i>Usted puede aplicar al equipo cualquiera de los siguientes pictogramas:</i>
<i>français</i>	<b><i>Symboles utilisés sur l'instrument</i></b> <i>Tous les symboles suivants peuvent être utilisés sur l'instrument :</i>
<i>italiano</i>	<b><i>Contrassegni utilizzati sullo strumento</i></b> <i>Qualsiasi dei seguenti contrassegni può essere utilizzato sullo strumento:</i>
<i>Nederlands</i>	<b><i>Markeringen op het Instrument</i></b> <i>De volgende markingen kunnen op het instrument gebruikt worden:</i>
<i>português</i>	<b><i>Marcações usadas no aparelho</i></b> <i>Qualquer uma das marcações seguintes podem ser usadas no aparelho:</i>
<i>suomea</i>	<b><i>Laitteessa käytetyt merkit</i></b> <i>Laitteessa voidaan käyttää seuraavia merkkejä:</i>
<i>svenska</i>	<b><i>Apparatens symboler</i></b> <i>Följande symboler kan finnas på apparaten:</i>

This symbol indicates that there is a potential hazard and that the user must look in the guide for an explanation of the hazard and how to avoid it.

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This symbol indicates the operational on/off switch.

---

<i>dansk</i>	Dette symbol angiver betjenings-tændt/sluk-kontakt.
<i>Deutsch</i>	Dieses Symbol kennzeichnet den Bereitschaftsschalter.
<i>español</i>	Este símbolo indica el interruptor operacional con./desc.
<i>français</i>	Ce symbole représente l'interrupteur d'arrêt/marche.
<i>italiano</i>	Questo simbolo indica l'interruttore acceso/spento opzionale.
<i>Nederlands</i>	Dit symbool geeft de aan/uit schakelaar aan.
<i>português</i>	Este símbolo indica o interruptor de ligar/desligar o funcionamento.
<i>suomea</i>	Tämä symboli tarkoittaa päälle/pois -käyttökytkintä.
<i>svenska</i>	Symbolen visar apparatens till/från-knapp.

---

This symbol indicates alternating current.

---

<i>dansk</i>	Dette symbol angiver vekselstrøm.
<i>Deutsch</i>	Dieses Symbol kennzeichnet Wechselstrom.
<i>español</i>	Este símbolo indica la existencia de corriente alterna.
<i>français</i>	Ce symbole représente le courant alternatif.
<i>italiano</i>	Questo simbolo indica corrente alternata.
<i>Nederlands</i>	Dit symbool geeft wisselstroom aan.
<i>português</i>	Este símbolo indica corrente alterna.
<i>suomea</i>	Tämä symboli merkitsee vaihtovirtaa.
<i>svenska</i>	Symbolen indikerar växelström.

This symbol indicates that there is a potential hazard due to a strong magnetic field and that persons wearing heart pacemakers or other metallic implants should remain at least 0.6 m away from the furnace.

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## Warning messages

This section contains translations of the warnings used in this guide.

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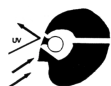
---

**Caution:**

The term CAUTION alerts you to situations that could result in **serious damage to the instrument** or other equipment.

Details about these circumstances are described in a message similar to this one.

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**Warning: Explosive Atmosphere**

This instrument is not designed for operation in an explosive atmosphere.

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**Warning: Electrical Hazard**

Any interruption of the protective conductor inside or outside the instrument or disconnection of the protective conductor (earth/ground) terminal is likely to make the instrument dangerous.

Intentional interruption is prohibited.

**Lethal voltages are present in the instrument**

- Even with the power switch off, line power voltages can still be present within the instrument.
- When the instrument is connected to line power, terminals may be live, and opening covers or removing parts (except those to which access can be gained without use of a tool) is likely to expose live parts.
- Capacitors inside the instrument may still be charged even if the instrument has been disconnected from all voltage sources.

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**A**

- AAAnalyst 800
  - concept 8-3
  - features 8-3
  - system components 8-4
  - technical data 8-13
- acetylene 2-15
  - connecting 3-7
- acetylides 2-18
- air
  - connecting 3-7
- air filters
  - changing 5-9
- argon
  - connecting 3-10
- automatic circuit breaker
  - resetting 5-7

**B**

- burner drain system
  - connecting 3-14
- burner gases
  - acetylene 2-15
  - connecting 3-6
- burner system
  - drain system 2-17
  - safe operation 2-16

**C**

- changing air filters 5-9
- changing fuses 5-7
- checklist
  - general maintenance 5-4
  - spectrometer maintenance 5-4
- circuit breaker 4-3
  - resetting 5-7
- cleaning windows 5-5
- compressed gases 2-10

## connecting

- acetylene 3-7
- air 3-7
- argon 3-10
- burner drain system 3-14
- burner gases 3-6
- furnaces gases 3-9
- nitrogen 3-10
- nitrous oxide 3-7
- special gas 3-12
- connections
  - electrical 3-18
- connectors on the spectrometer
  - electrical 8-10
  - gas and cooling 8-8
- cooling system 7-15
  - connecting and disconnecting
    - water hoses 3-15
  - filling 3-17
  - furnace cooling 7-15
  - installing 3-15
  - spectrometer cooling 7-15
- correct use of the instrument 1-7, 2-3
- customer service
  - catalog service 1-3
  - company name and address 1-3
  - supplies, replacement parts, and accessories 1-3
- cyanide solutions 2-18

**D**

- deuterium lamp
  - replacing 5-10
- dimensions of the system 7-4
- disposing of
  - contents of waste containers 2-7
  - defective lamps 2-7
- drain system 2-17

**E**

- electrical components
  - maintaining 5-7
- electrical connections 3-18
- electrical protection 1-4, 8-15
- electrical requirements 7-8
- electrical safety 1-8
- electrodeless discharge lamps
  - installing and connecting 4-8
- electromagnetic compatibility 1-5
- electrothermal atomization
  - hazards 2-20
  - safety practices 2-19
- emergency shutdown 4-14
- environment
  - operating conditions 1-10
  - storage conditions 1-10
- environmental conditions 7-9

**F**

- FIAS and mercury/hydride systems
  - hazards 2-22
  - safety practices 2-21
- filling the cooling system 3-17
- flame
  - safe operation 2-16
- flame atomization
  - safe operation of the flame 2-16
  - safe use of burner gases 2-13
  - safety practices 2-13
- flashbacks 2-17
- fume ventilation 7-10
- furnace autosampler
  - installing 3-14
- furnace gases
  - connecting 3-9

fuses  
changing 5-7

## G

gas hazards 2-10  
gas supply requirements 7-11  
air 7-13  
burner gases 7-12  
furnace gases 7-14

## H

hazards with electrothermal atomization  
high temperatures 2-20  
toxic products 2-20  
UV radiation 2-20  
hazards with FIAS and mercury/hydride systems  
hazardous chemicals 2-22  
high temperatures 2-22  
toxic products 2-22  
hazards with flame atomization  
acetylides 2-18  
cyanide solutions 2-18  
high temperatures 2-18  
toxic combustion products 2-18  
UV radiation 2-18  
heart pacemakers 2-8  
high temperatures 2-9  
hollow cathode lamps  
installing 4-8  
hygiene 2-4

## I

IEC 1010 compliance 1-4  
installation procedures 3-3  
installing  
cooling system 3-15  
lamps 4-7  
sample trays 4-11

installing and connecting lamps 4-8  
installing furnace autosampler 3-14

## L

laboratory  
hygiene 2-4  
ventilation 2-4  
laboratory requirements  
cooling system 7-15  
environmental conditions 7-9  
fume ventilation 7-10  
gas supplies 7-11  
lamp compartment 4-7  
lamps  
installing 4-7  
installing and connecting 4-8  
removing 4-10

## M

magnetic field 2-8  
safe radius 7-7  
maintaining electrical components 5-7  
maintenance checklists 5-4  
maintenance procedures 5-5  
markings on the instrument 1-12  
moving the spectrometer 3-4

## N

nitrogen  
connecting 3-10  
nitrous oxide  
connecting 3-7

## O

operating conditions 1-10  
operational on/off switches 8-12  
optical system 8-6

## P

parts and supplies 6-1  
parts provided 6-3  
Perkin-Elmer service 5-3

## R

Reconfigure utility 4-4  
references for laboratory safety practice 2-23  
regulatory information  
electromagnetic compatibility 1-5  
removing lamps 4-10  
replacement parts 6-5  
catalog service 6-5  
replacing air filters 5-9  
replacing fuses 5-7  
replacing the deuterium lamp 5-10  
resetting the circuit breaker 5-7

## S

safe handling of chemicals 2-5  
sodium tetrahydroborate 2-6  
safe use of burner gases  
acetylene 2-15  
safety  
compressed gases 2-10  
correct use of the instrument 1-7, 2-3  
cyanide solutions 2-18  
electrical safety 1-8  
handling chemicals 2-5  
high temperatures 2-9  
magnetic field 2-8  
safe radius 7-7  
operation of the flame 2-16  
UV radiation 2-8  
safety conventions 1-8  
see also symbols

safety practices  
  electrothermal atomization  
    2-19  
  FIAS and mercury/hydride  
    systems 2-21  
  flame atomization 2-13  
  references 2-23  
sample trays  
  installing 4-11  
shutting down the system 4-12  
sodium tetrahydroborate 2-6  
space requirements 7-4  
special gas  
  connecting 3-12  
spectrometer  
  moving 3-4  
spectrometer maintenance 5-5  
storage conditions 1-10  
switching on the system 4-3  
symbols  
  used in the handbook 1-6  
  used on the instrument 1-11  
  translations T-3  
system  
  shutting down 4-12

## T

technical data  
  spectrometer 8-13  
translations  
  symbols used on the instru-  
    ment T-3  
  warning messages T-11  
troubleshooting 5-3

## U

UV radiation 2-8

## V

ventilation, laboratory 2-4

## W

warning markings on the instru-  
  ment 1-12  
waste disposal  
  contents of waste containers  
    2-7  
  disposing of defective lamps  
    2-7  
windows  
  cleaning 5-5

## Z

Zeeman magnetic field 2-8, 7-7