Agilent 7100 Capillary Electrophoresis System

User Manual

Agilent Technologies

Notices

© Agilent Technologies, Inc. 2009 - 2019

No part of this manual may be reproduced in any form or by any means (including electronic storage and retrieval or translation into a foreign language) without prior agreement and written consent from Agilent Technologies, Inc. as governed by United States and international copyright laws.

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Research Use Only

Warranty

The material contained in this document is provided “as is,” and is subject to being changed, without notice, in future editions. Further, to the maximum extent permitted by applicable law, Agilent disclaims all warranties, either express or implied, with regard to this manual and any information contained herein, including but not limited to the implied warranties of merchantability and fitness for a particular purpose. Agilent shall not be liable for errors or for incidental or consequential damages in connection with the furnishing, use, or performance of this document or of any information contained herein. Should Agilent and the user have a separate written agreement with warranty terms covering the material in this document that conflict with these terms, the warranty terms in the separate agreement shall control.

Technology Licenses

The hardware and/or software described in this document are furnished under a license and may be used or copied only in accordance with the terms of such license.

Restricted Rights Legend

If software is for use in the performance of a U.S. Government prime contract or subcontract, Software is delivered and licensed as “Commercial computer software” as defined in DFAR 252.227-7014 (June 1995), or as a “commercial item” as defined in FAR 2.101(a) or as “Restricted computer software” as defined in FAR 52.227-19 (June 1987) or any equivalent agency regulation or contract clause. Use, duplication or disclosure of Software is subject to Agilent Technologies’ standard commercial license terms, and non-DOD Departments and Agencies of the U.S. Government will receive no greater than Restricted Rights as

defined in FAR 52.227-19(c)(1-2) (June 1987). U.S. Government users will receive no greater than Limited Rights as defined in FAR 52.227-14 (June 1987) or DFAR 252.227-7015 (b)(2) (November 1995), as applicable in any technical data.

Safety Notices

CAUTION

A **CAUTION** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a **CAUTION** notice until the indicated conditions are fully understood and met.

WARNING

A **WARNING** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a **WARNING** notice until the indicated conditions are fully understood and met.

In This Guide...

This guide contains information to install and operate the Agilent 7100 Capillary Electrophoresis System.

1 Introduction

This chapter gives an introduction to the instrument, instrument overview and internal connectors.

2 Site Requirements and Specifications

This chapter provides information on environmental requirements, physical and performance specifications.

3 Installing the Instrument

In this chapter you will find information about how to unpack and to set up the Agilent CE and information about the configuration of Instrument.

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6 Troubleshooting and Diagnostics

This chapter gives an overview about the troubleshooting and diagnostic features and the different user interfaces.

7 Hardware Information

This chapter describes the instrument in more detail on hardware and electronics.

8 Maintenance

This chapter describes the maintenance of the CE instrument.

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This chapter provides information on parts for maintenance.

10 Appendix

This chapter provides additional information on safety, legal and web content associated with your 7100 CE instrument.

Contents

1	Introduction	13
	Introduction to the Agilent 7100 Capillary Electrophoresis System	14
	Getting Help	14
	Where to Find Information	15
	What Learning Products are Available?	16
2	Site Requirements and Specifications	17
	Site Requirements	18
	Power Consideration	18
	Power Cords	19
	Bench Space	20
	Environment	20
	Physical Specifications	21
	Performance Specifications	22
	Platzbedarf	25
	Stromversorgung	25
	Netzkabel	26
	Platzbedarf	27
	Umgebungsbedingungen	27
	Abmessungen und Gewicht	28
	Leistungsspezifikationen	29
	Exigences d'installation	32
	Remarques à propos de l'alimentation	32
	Câbles d'alimentation	33
	Encombrement	34
	Environnement	34
	Caractéristiques physiques	35
	Caractéristiques de performance	36

Contents

Requisiti ambientali	39
Alimentazione	39
Cavi di alimentazione	40
Spazio necessario	41
Ambiente	41
Specifiche fisiche	42
Specifiche delle prestazioni	43
Requisitos de las instalaciones	46
Consideraciones sobre la corriente	46
Cables de alimentación	47
Espacio en el banco	48
Entorno	48
Especificaciones físicas	49
Especificaciones de rendimiento	50
設置について	53
電源について	53
電源コード	54
作業台スペース	55
環境条件	55
物理的仕様	56
性能仕様	57
场地要求	60
电源要求	60
电源线	61
工作台位置	62
环境	62
物理规格	63
性能规格	64
3 Installing the Instrument	67
Unpacking the Agilent CE Instrument	68
Damaged Packaging	68
Unpacking the CE Instrument	68

Delivery Checklist	69
Installing the Instrument	71
Physical Connection	71
4 LAN Configuration	75
What You Have to Do First	76
TCP/IP Parameter Configuration	77
Configuration Switch	78
Initialization Mode Selection	79
Link Configuration Selection	83
Storing the Settings Permanently with Bootp	84
Manual Configuration	85
With Telnet	86
Agilent Bootp Service	89
Location of the Media Access Control (MAC) Address	90
5 Using the Agilent 7100 Capillary Electrophoresis System	91
System Definition	92
Controlling the Agilent CE Instrument	93
Before You Start the Agilent CE Instrument	93
Starting Up the Agilent CE Instrument	93
Starting Your Computer and the Agilent ChemStation Software	95
Starting the Agilent ChemStation Software the First Time	96
Working with the CE Diagram	98
Common Software Toolbar	98
Software Toolbar	99
Graphical User Interface Explanation	100
Menus in the CE and DAD Diagram	102
System Vialtable	104
How to Prepare your Agilent CE Instrument for Analysis	105
Capillary Electrophoresis, CE	106
How to Use Your Agilent CE Instrument for Capillary Electrophoresis Separations	106

Contents


Agilent CE Method Parameters	106
Vials	110
Cassette Temperature	111
High Voltage System	112
Posttime and Stoptime	114
Replenishment	114
Preconditioning	115
Injection	117
Timetable	119
Detector Parameters	121
Switching the Lamp On	124
Preparing the Replenishment System if Needed	125
Preparing Vials for Buffer and Sample	129
Installing a Capillary in the Alignment Interface	137
Installing a Capillary in an Empty Capillary Cassette	141
Inserting the Cassette	145
Fraction Collection	148
Capillary Electrophoresis, Plus High Pressure CE+p	157
Requirements	157
Outlet Lift Connection to the Tubing System	159
Preparing the Agilent CE Instrument for High Pressure Use	161
Operation and Function in CE+p Mode	164
Injection Options with CE+p	167
Timetable Options with CE+p	167
CE+p Control Using the CE Diagram	168
Capillary Electrochromatography, CEC	169
How to Use Your Agilent CE Instrument for Capillary Electrochromatography	169
Preparing the Agilent CE Instrument for High Pressure Use	171
CEC-Specific GUI Explanation	173
Injection Options with CEC	176
Timetable Options with CEC	177
CEC Control Using the CE Diagram	178

Running the CEC Analysis	179
6 Troubleshooting and Diagnostics	181
CE Problem Solving	182
Power On Fail	182
Leaks	183
Blocked Replenishment Needle	185
Leakage Current	185
Broken Capillary	186
Contaminated Insulation Plate	186
Empty Injection Vial	187
Dirt in the Vials	187
Air Bubble in Injection Vial or Run Buffer Vials	187
Air Bubble in Capillary	187
Problems with the Replenishment System and Buffers Containing SDS	188
Agilent Lab Advisor Software	189
Configuration	189
7 Hardware Information	191
Identifying Instrument Components	192
7100 Capillary Electrophoresis Instrument Electronics	193
Firmware	196
Resident System	196
Main System	196
Firmware Updates	197
Diode Array Detector	199
DAD Control	200
Digital to Analog Converter	201
Changing the Fuses for the Power Supplies	202
Installing Drainage Tubing	203
External Water Bath for Tray Cooling (Optional)	204
Analog to Digital Converter	206

8	Maintenance	209
	Standard Operating Procedures for Maintenance of Your Agilent CE Instrument	210
	Overview of Maintenance	211
	Early Maintenance Feedback (EMF)	214
	EMF Counters	214
	Using the EMF Counters	214
	Setting the EMF Limits	215
	Cleaning the Electrodes, Pre-punchers and Insulation Plate	216
	Prepare the Agilent CE Instrument	217
	Remove the Covers	218
	Accessing the Electrodes	219
	Accessing the Liquid Handling Module	221
	Accessing and Removing the Pre-punchers	223
	Cleaning the Electrodes	225
	Cleaning the Pre-punchers	226
	Cleaning the Insulation Plate	227
	Reinstalling the Pre-punchers	228
	Reinstalling the Electrodes and the Insulation Plate	230
	Reinstall the Capillary Cassette	235
	Reinstall the Front Cover	235
	Maintenance of the Replenishment System	236
	Changing the Buffer Composition in the Replenishment System	237
	The Replenishment System Will Not Be Used for Some Time	238
	Further Action	239
	Operations for Replenishment Maintenance	239
	Changing the Air Inlet Filter	242
	Accessing the Air Filter	242
	Changing the Lamp	244
	Removing the Existing Lamp	245
	Inserting a New Lamp	247
	Further Action	247
	Cleaning the Instrument	248

9	Parts and Materials for Maintenance	249
	Overview of Maintenance Parts	250
10	Appendix	251
	General Safety Information	252
	Operation	252
	Allgemeine Sicherheitsinformation	255
	Informations générales de sécurité	258
	Informazioni generali sulla sicurezza	261
	Información de seguridad	264
	安全に関する一般的な情報	267
	一般安全信息	270
	Setting Up a Test Sample Run	272
	Method Parameter CE (Default Settings)	273
	Method Parameter DAD	275
	Chemical and Biological Safety	276
	The Waste Electrical and Electronic Equipment (WEEE) Directive (2002/96/EC)	277
	Radio Interference	278
	Sound Emission	279
	Legal Notice	280
	Agilent Technologies on Internet	281
	Index	283

Contents

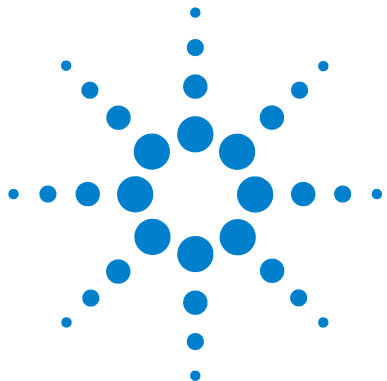1 Introduction

Introduction to the Agilent 7100 Capillary Electrophoresis System	14
Where to Find Information	15
What Learning Products are Available?	16

This chapter gives an introduction to the Agilent 7100 Capillary Electrophoresis System and where to find specific information or help.

Introduction to the Agilent 7100 Capillary Electrophoresis System

This handbook gives an overview of your Agilent 7100 Capillary Electrophoresis System. It is designed to get you started working with the system. We strongly advise you to make extensive use of the online help when working with the system. There you find detailed reference and task information which complements the overview given in this guide. Refer to “[Getting Help](#)” later in this section for information on using the online help of the ChemStation. Refer to “[Where to Find Information](#)” for help in finding information on certain tasks.

Getting Help

You can get help by choosing the **Help** item from the top menu. This puts you into the help index. From there you can browse through the information by selecting the appropriate jumps (underlined words). To access specific information in the help:

When you are in the Agilent ChemStation or Lab Advisor software:

- select a context-sensitive element and press F1.

When you are in the online help:

- search for a specific keyword using the search command,
- select a jump - a jump is an underlined word you can select to go to a topic related to the word, or
- select a word or phrase with a dotted underline to see its definition. The online help also contains online information on how to use help.

Where to Find Information

Table 1 Where to find information

Task	Refer to
• Installation of the Agilent CE System, installing the external water bath, installing the high sensitivity cell	"The Core Agilent CE Instrument" " Installing the Instrument " on page 67
• Setting up the Agilent CE System for CE analysis, CE+p	Chapter " Using the Agilent 7100 Capillary Electrophoresis System " on page 91 "
• How to use the Agilent CE System in the CE+p mode	" Capillary Electrophoresis, Plus High Pressure CE+p " on page 157
• How to use the Agilent CE System for capillary electrochromatography	" Capillary Electrochromatography, CEC " on page 169
• How to maintain the Agilent CE System	Chapter " Maintenance " on page 209
• Setting up a sequence	Online help " Understanding Your ChemStation "
• Data analysis	Online help " Understanding Your ChemStation "
• Setting up a report	Online help " Understanding Your ChemStation "
• Spectral library search	" ChemStation Applications " handbook and online help
• Setting up customized reports	" ChemStation Applications " handbook
• Safety information	Safety information in " Appendix " on page 251
• Learning ChemStation concepts	" Understanding Your ChemStation " handbook
• Learning the theory of the technique	<i>High Performance Capillary Electrophoresis: An Introduction</i> The CE Partner CD-ROM: " <i>The comprehensive, interactive tool for beginners and advanced users of CE</i> "
• Application	Agilent web page (http://www.agilent.com), search for Application notes


What Learning Products are Available?

The Agilent CE System is supplied with the following:

- User's Guide
- High Performance Capillary Electrophoresis: An Introduction

The Agilent ChemStation and Lab Advisor/Instrument Utilities software DVD is supplied with the following electronic documents:

- Installing Your Agilent ChemStation for LC and CE Systems
- Installing and Understanding Your Spectra Module
- Understanding Your Agilent ChemStation
- Online help
- Understanding Your Agilent ChemStation for CE Systems

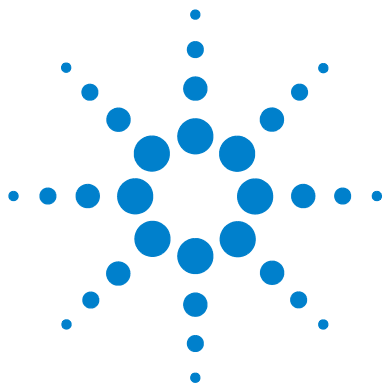2 Site Requirements and Specifications

Site Requirements	18
Physical Specifications	21
Performance Specifications	22
Platzbedarf	25
Abmessungen und Gewicht	28
Leistungsspezifikationen	29
Exigences d'installation	32
Caractéristiques physiques	35
Caractéristiques de performance	36
Requisiti ambientali	39
Specifiche fisiche	42
Specifiche delle prestazioni	43
Requisitos de las instalaciones	46
Especificaciones físicas	49
Especificaciones de rendimiento	50
設置について	53
物理的仕様	56
性能仕様	57
场地要求	60
物理规格	63
性能规格	64

This chapter provides information on the environmental requirements, as well as physical and performance specifications.

Site Requirements

A suitable environment is important to ensure optimal performance of the 7100 Capillary Electrophoresis instrument.

Before you begin installation, check carefully that the place you have chosen meets the requirements below.

Power Consideration

The instrument power supply has wide ranging capabilities and accepts any standard line voltage in the range mentioned in [Table 2](#) on page 21. Consequently, there is no voltage selector on the back of the detector.

WARNING

The module is partially energized when switched off, as long as the power cord is plugged in.

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- Make sure that it is always possible to access the power plug.
- Remove the power cable from the instrument before opening the cover.
- Do not connect the power cable to the instrument while the electronic box in the back of the instrument is opened or side panels are removed.

WARNING

Incorrect line voltage on the instrument

Shock hazard or damage to your instrument can result if the device is connected to a line voltage that is higher than specified.

- Connect your instrument to the specified line voltage.
-

CAUTION

Inaccessible power plug.

In the case of an emergency it must be possible to disconnect the instrument from the power line at any time.

- Make sure the power connector of the instrument can be easily reached and unplugged.
 - Provide sufficient space behind the power socket of the instrument to unplug the cable.
-

Power Cords

Different power cords are delivered with the instrument. The female end of all power cords is identical. It plugs into the power-input socket at the rear. The male end of each power cord is different and designed to match the wall socket of a particular country or region.

WARNING

Absence of a ground connection or use of an unspecified power cord

The absence of a ground connection or the use of an unspecified power cord can lead to an electric shock or short circuit.

- Never operate your instrument from a power outlet without a ground connection.
 - Never use a power cord other than the Agilent Technologies power cord designed for your region.
-

WARNING

Use of unsupplied cables

Using cables not supplied by Agilent Technologies can lead to damage of the electronic components or personal injury.

- Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.
-

Bench Space

The instrument dimensions and weight mean that the instrument can be placed on almost any desk or laboratory bench. It needs an additional 2.5 cm (1.0 inch) of space on either side and approximately 8 cm (3.1 inches) in the rear for air circulation and electric connections.

If the Agilent 7100 Capillary Electrophoresis instrument is to be placed on a bench, make sure that the bench is designed to bear the weight of the system.

The instrument must be operated in a upright position.

Environment

Your instrument will work within the specifications for ambient temperatures and relative humidity described in [Table 2](#) on page 21.

CAUTION

Condensation within the module

Condensation will damage the system electronics.

- Do not store, ship or use your module under conditions where temperature fluctuations could cause condensation within the module.
 - If your module was shipped in cold weather, leave it in its box and allow it to warm slowly to room temperature to avoid condensation.
-

Physical Specifications

Table 2 Physical specifications

Type	Specification	Comments
Weight	35 kg (77.2 lbs)	
Dimensions (width × depth × height)	350 x 510 x 590 mm (13.8 x 20.1 x 23.2 inches)	
Line voltage	100–240 VAC ± 10%	Wide-ranging capability
Line frequency	50 or 60 Hz ± 5%	
Power consumption	350 VA / 300 W / 1024 BTU/h	Maximum
Ambient operating temperature	5–40 °C (41–104 °F)	
Ambient non-operating temperature	-40 to 70°C (-40 to 158 °F)	
Humidity	below 80% at 31 °C (87.8 °F)	Non-condensing
Operating altitude	Up to 2000 m (6500 ft)	
External cooling	max. 0.5 bar (7.2 psi), max. 50 °C (122 °F)	Waterbath
External pressure	2-12 bar (29-174 psi)	Oil-free air or nitrogen
Safety standards: IEC, CSA, UL	Installation category II, Pollution degree 2	For indoor use only
Housing	All materials are recyclable	

For applications run in CEC mode (Capillary Electrochromatography) or CE+p mode (Capillary Electrophoresis mode with optional usage of higher pressure), an external pressure supply of oil-free air or nitrogen with a maximum pressure of 15 bar (218 psi) can be used. A respective male adapter that fits into the instrument's female adapter and PTFE tubing are part of the shipment.

NOTE

The Agilent 7100 Capillary Electrophoresis System is designed to operate in a typical electromagnetic environment where RF transmitters, such as mobile phones, should not be used in close proximity.

Performance Specifications

Table 3 Performance specifications for the Agilent 7100 Capillary Electrophoresis System

Type	Specification	Comments
Safety features	Extensive diagnostics, error detection and display, current leak detection; low current limit Liquid leak sensor; safety sensors on door and cover disabling high voltage; vial sensor;	
Control and data evaluation	Agilent control software (e.g. ChemStation, OpenLAB CDS ChemStation Edition, MassHunter)	
Communication	LAN, controller-area network (CAN), RS-232C, APG remote: ready, start, stop and shut-down signals. USB (future use). Analog out and in.	0 to 1 V
GLP features	Early maintenance feedback (EMF), electronic records of maintenance and errors	
Diagnostic functions	RFID tag for lamp, early maintenance feedback, supported by Agilent Lab Advisor software with integrated diagnosis test suite	
Pressure system	<ul style="list-style-type: none"> • Programmable with -100 to +100 mbar on inlet • Flushing with 1 bar or with high pressure 2–12 bar • Vial pressurization with high pressure 2–12 bar on inlet and/or outlet 	
Injection modes	<ul style="list-style-type: none"> • Self correcting injection system with injection from inlet or outlet • Programmable range • Pressure • Electrokinetic 	up to 10,000 seconds –100 to +100 mbar –30 to +30 kV
Replenishment	<ul style="list-style-type: none"> • Satellite station for refilling of inlet or outlet buffer vials with fresh buffer for automatic, continuous operation. • Selectable buffer levelling. 	
Vials	<ul style="list-style-type: none"> • 100 µL (polypropylene or glas) with resealing snap caps • 1 mL (polypropylene) with resealing snap caps • 2 mL (glas) with resealing snap caps 	

Table 3 Performance specifications for the Agilent 7100 Capillary Electrophoresis System

Type	Specification	Comments
Raw data channels	Detector signals and spectrum, voltage, current, leak current, power, cassette temperature, pressure, lamp voltage and detector temperature	
Electrophoresis power	<ul style="list-style-type: none"> • Voltage range • Current • Power • Operation under constant voltage, current or power • Programmable polarity switch 	setable 0 to ± 30 kV supply setable 0–300 μ A setable 0–6 W
CE specific software	<ul style="list-style-type: none"> • Mobility report, time corrected areas, pI calibration and bio polymer size calibration 	

Table 4 Performance specifications for the Agilent 7100 Capillary Electrophoresis System, Autosampler/ fraction collector

Type	Specification	Comments
Autosampler/ fraction collector	50-position carousel All vials are randomly accessible from inlet or outlet end of capillary Temperature control with external waterbath with vial temperature from 10–40 °C (non-condensing conditions, minimum waterbath temperature +1 °C)	

Table 5 Performance specifications for the Agilent 7100 Capillary Electrophoresis System, Detector

Type	Specification	Comments
Detector	Real time UV-Visible diode-array detector (190–600 nm)	
	Temperature controlled	
Wavelength accuracy	1 nm	
Response time	0.063 to 16 s	
Light source	Prealigned deuterium lamp	

2 Site Requirements and Specifications

Performance Specifications

Table 5 Performance specifications for the Agilent 7100 Capillary Electrophoresis System, Detector

Type	Specification	Comments
Signals	Up to eight signals simultaneously, full spectral acquisition with Agilent ChemStation	
Sensitivity	1 μ M 4-hydroxy-aceto-phenon injected at 50 mbar*5 sec, 3 x 50 μ m bubble cell capillary, signal/noise >6* (50 mM borate buffer pH 9.3, 25 kV)	
Baseline noise	<50 μ AU (at 2 sec response time)*	
Linear dynamic range	1x10 ⁴ (3x50 μ m bubble cell capillary)*	

* typical value

Table 6 Performance specifications for the Agilent 7100 Capillary Electrophoresis System, Capillary Cassette

Type	Specification	Comments
Capillary cassette	High-speed forced-air cooler with Peltier element	
Temperature range	10 °C below ambient, up to 60 °C (min. 10 °C cassette temperature)	
Minimum total capillary length	31 cm	
Capillary compatibility	365 μ m o.d.	

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3 Installing the Instrument

Unpacking the Agilent CE Instrument 68
Installing the Instrument 71

In this chapter you will find information on

- how to unpack the Agilent CE instrument
- how to set up the Agilent CE instrument

Information on how to install the Agilent ChemStation software or on how to install the Lab Advisor/Instrument Utilities software can be found in separate documentation located on the respective software DVD.

Unpacking the Agilent CE Instrument

Damaged Packaging

If the delivery packaging shows signs of external damage, please call your Agilent Technologies sales and service office immediately. Inform your service representative that the instrument may have been damaged during shipment.

CAUTION

"Defective on arrival" problems

If there are signs of damage, please do not attempt to install the module. Inspection by Agilent is required to evaluate if the instrument is in good condition or damaged.

- Notify your Agilent sales and service office about the damage.
 - An Agilent service representative will inspect the instrument at your site and initiate appropriate actions.
-

Unpacking the CE Instrument

The Agilent CE instrument is heavy and should be carried by two people.

- 1** Use the recessed handles on both sides to carry the Agilent CE instrument.
- 2** Open the box in which the Agilent CE instrument was shipped.
- 3** Pull the Agilent CE instrument out of the box and position it in a suitable location.

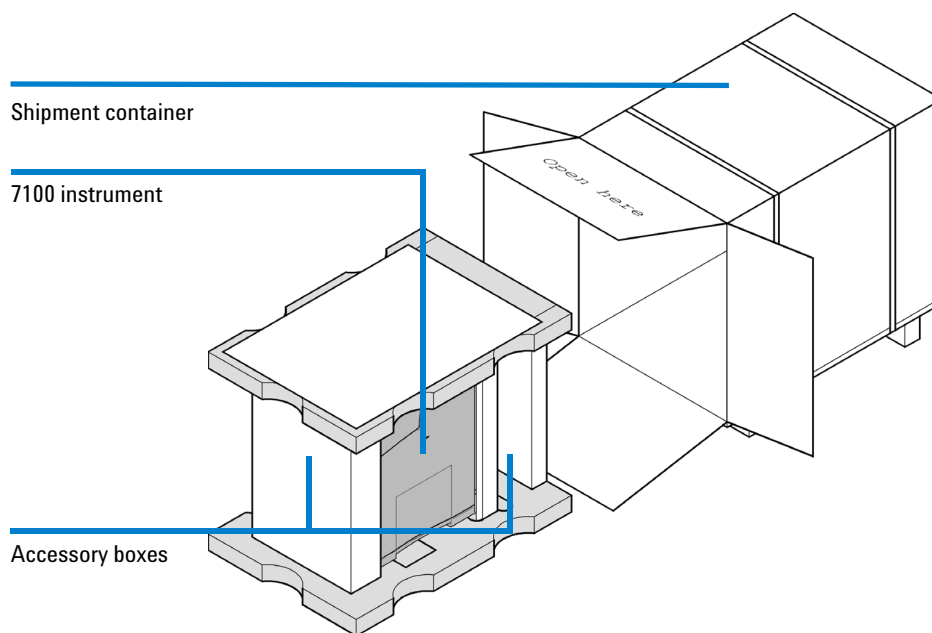
Figure 1 Unpacking the Agilent CE instrument

Delivery Checklist

Ensure all parts and materials have been delivered with the instrument. Please report missing or damaged parts to your local Agilent Technologies sales and service office.

Accessories

The accessories shipped with the Agilent CE instrument are in additional boxes on right and left in the instrument box. Check that you have received everything using [Table 6](#).

3 Installing the Instrument

Unpacking the Agilent CE Instrument

Table 6 Accessories

Description	Part number	Quantity
Accessory kit	G7100-68705	1
– HPCE startup capillary kit	1600-68700	1
– Detector interface, green, 50 µm	G7100-60210	1
– Detector interface, red, extended 50 µm	G7100-60230	1
– Coupler	0100-1633	2
– O-ring 25.12-mm-id 1.78-mm-XSECT-DIA	0905-1163	1
– O-ring 3-mm	0905-1196*	5
– Container	1540-0307	1
– Fuse T6.3AH, 250 VAC	2110-1018	2
– Fuse T6.3AL, 250 VAC	2110-0623	2
– Fuse T3.15AL, 250 VAC	2110-1417	4
– Air filter	3150-0619	1
– Tubing assembly	5063-6527	1
– Screwdriver, slot shaft	8710-2607	1
– Screwdriver, Pozidriv shaft	8710-0899	1
– 12 mm socket wrench	8710-2076	1
– Lift tubing assy	G1600-67003	1
– Vial crimp/snap polypropylene 1 mL 100/PK	5182-0567	1
– Snap caps, polyurethane 100/PK	5181-1512	1
– Cable, remote	5061-3378	1
– Bottle, 500 mL, electrolyte, pressure plus	9301-6531	2
– Bottle, 100 mL, electrolyte, pressure plus	9301-6532	1
– CE System User’s Guide	G7100-90000	1
– Glass filter, solvent inlet	5041-2168	1
– Electrode alignment assy	G7100-60107	1
– LAN cable (crossover)	5023-0203	1
CE Installation Qualification Kit	5063-6514	1

*Part number 0905-1196 is not orderable. Please re-order part number 5062-8544 (5/PK)

Installing the Instrument

Physical Connection

WARNING

Module is partially energized when switched off, as long as the power cord is plugged in.

Repair or maintenance work on the module can lead to personal injuries, for example a shock hazard, when the cover is opened and the module is connected to the power.

- Ensure that line power cord is disconnected before setting up your Agilent CE instrument.

In general you have to go through the following steps.

- 1 Unpack the Agilent CE instrument and position it in a suitable location on your bench, see [“Unpacking the Agilent CE Instrument”](#) on page 68.
- 2 Connect the cables at the rear of the Agilent CE instrument.
- 3 Install the bottles in the replenishment system, see [“Preparing the Replenishment System if Needed”](#) on page 125.

Connecting Cables

The Agilent CE instrument has the following connectors on the rear.

- All electrical connectors, such as line power, LAN connector or fuses, are located on the right-hand side of the electronic box on the back of the instrument.
- All connectors for liquids and external pressure supply are located on the left-hand side of the electronic box.

- 1** Ensure that the line power switch on the front of the instrument is OFF.
- 2** Connect the following cables as shown in [Table 2](#) and in [Figure 2](#).
 - a** One power cable
 - b** LAN cable, which connects the Agilent CE instrument directly to the computer with the Agilent ChemStation installed or to the local network.

For a description of remaining items refer to the Hardware section.

- For more information on cable connection to 3rd party detectors or analog out, please refer to “[Hardware Information](#)” on page 191.

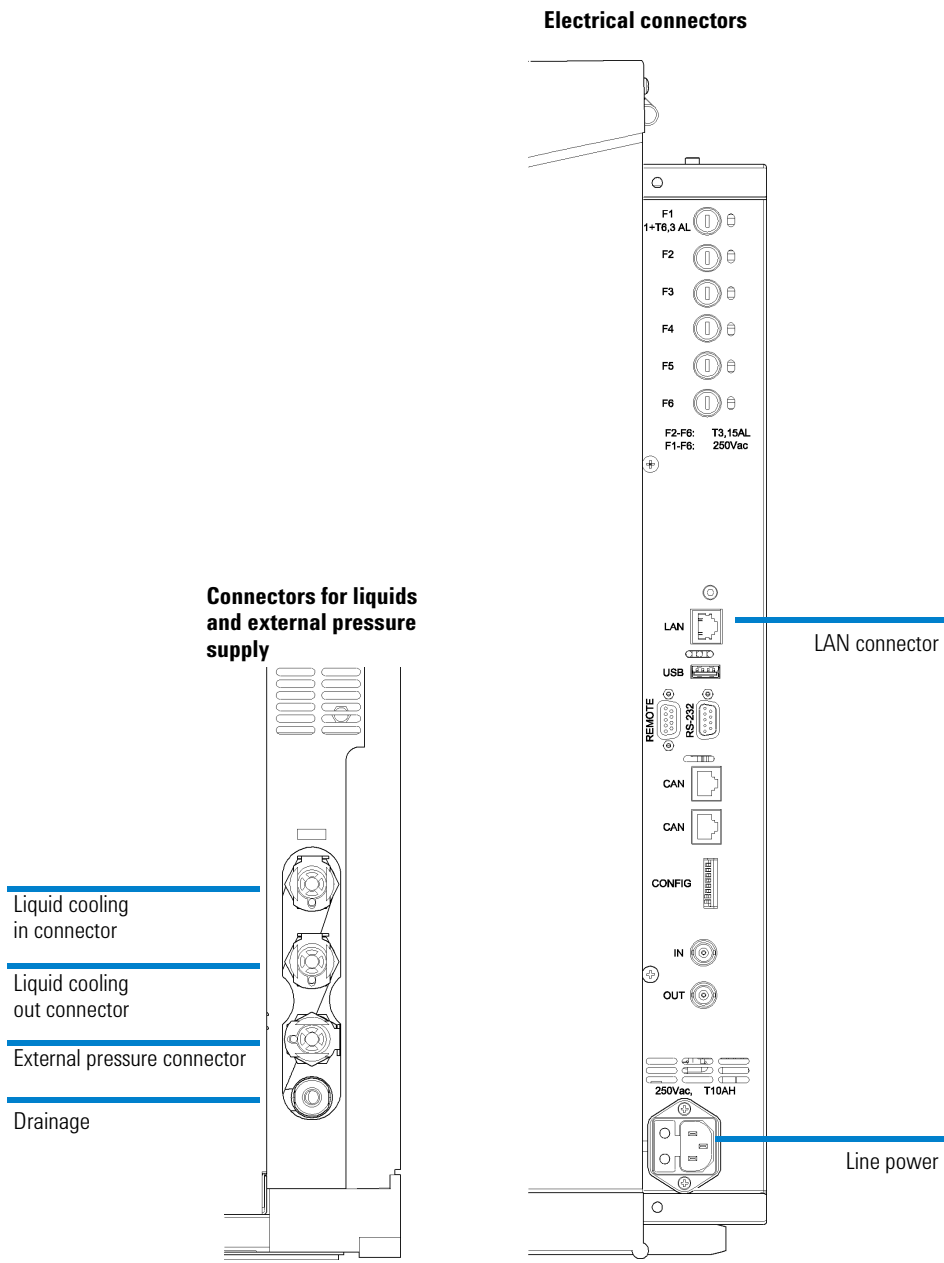
Figure 2 Connections on the rear of the instrument

3 Installing the Instrument

Installing the Instrument

4 LAN Configuration

What You Have to Do First	76
TCP/IP Parameter Configuration	77
Configuration Switch	78
Initialization Mode Selection	79
Link Configuration Selection	83
Storing the Settings Permanently with Bootp	84
Manual Configuration	85
Agilent Bootp Service	89
Location of the Media Access Control (MAC) Address	90

This chapter provides information on connecting the CE instrument to the Agilent ChemStation PC.

What You Have to Do First

The instrument has an on-board LAN communication interface. Please review the following options:

- 1 The instrument will be connected to the LAN: Ask your administrator to provide you with an IP address, a subnet mask address and a gateway address. Proceed with manual configuration (see [“Manual Configuration”](#) on page 85). You will have to initially establish a LAN connection using the default settings of the instrument and a crossover LAN cable (see [Table 9](#) on page 82).
- 2 The instrument and PC will be set up locally via a crossover cable or a hub: No addresses are needed. Proceed with establishing a LAN connection using the default settings of the instrument (see [Table 9](#) on page 82).
- 3 Identification of the instrument in your LAN by MAC address: If you plan to let a Bootp service (see [“Agilent Bootp Service”](#) on page 89) assign an IP address please continue with locating the Media Access (MAC) address and proceed with the respective initialization modes in [“Initialization Mode Selection”](#) on page 79 and the installation instruction in the "Installing your ChemStation for LC and CE" manual. Ask your administrator to provide you with an IP address, a subnet mask address and a gateway address.

In addition to LAN communication, the instrument allows serial communication access via the RS 232 cable to the COM port of the PC. However, the data transfer rate would be too low for running the instrument at full performance. This alternative data connection is only meaningful if the Lab Advisor software is for example used for troubleshooting reasons (see [“Agilent Lab Advisor Software”](#) on page 189).

Please see [“Starting Your Computer and the Agilent ChemStation Software”](#) on page 95 for information on configuration of the connection settings of the ChemStation software.

TCP/IP Parameter Configuration

To operate properly in a network environment, the LAN interface must be configured with valid TCP/IP network parameters. These parameters are:

- IP address
- Subnet mask
- Default gateway

The TCP/IP parameters can be configured by the following methods:

- by automatically requesting the parameters from a network-based BOOTP service (using the so-called bootstrap protocol), see [“Agilent Bootp Service”](#) on page 89
- by manually setting the parameters using Telnet

The LAN interface differentiates between several initialization modes. The initialization mode (abbreviation: 'init mode') defines how to determine the active TCP/IP parameters after power-on. The parameters may be derived from a Bootp cycle, non-volatile memory or initialized with known default values. The initialization mode is selected by the configuration switch, see [Table 7](#) on page 78.

Configuration Switch

The configuration switch can be accessed on the right rear side of the instrument. The 7100 CE is shipped with switches 7 and 8 set to ON. This means that the instrument is set to a fixed default IP: 192.168.254.11

NOTE

To perform any LAN configuration, SW1 and SW2 must be set to OFF.

Table 7 Factory default settings

Initialization ('init') mode	Using default, switches 7 and 8 set to ON.
Link configuration	Speed and duplex mode determined by auto-negotiation

Initialization Mode Selection

The following initialization (init) modes are selectable:

Table 8 Initialization mode switches

	SW6	SW7	SW8	Init mode
	OFF	OFF	OFF	Bootp
	OFF	OFF	ON	Bootp & Store
	OFF	ON	OFF	Using Stored
	OFF	ON	ON	Using Default

Bootp

When the initialization mode "Bootp" is selected, the 7100 CE tries to download the parameters from a Bootp server. The parameters obtained become the active parameters immediately. They are not stored in the non-volatile memory of the 7100 CE. Therefore, the parameters are lost with the next power cycle of the 7100 CE.

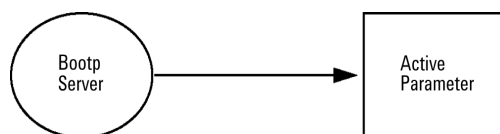
Figure 3 Bootp (principle)

Bootp & Store

When "Bootp & Store" is selected, the parameters obtained from a Bootp server become the active parameters immediately. In addition, they are stored in the non-volatile memory of the 7100 CE. Thus, after a power cycle they are still available. This enables a kind of "Bootp once" configuration of the 7100 CE.

Example: The user may not want to have a Bootp server be active in his network all the time. But on the other hand, he may not have any other configuration method than Bootp. In this case he starts the Bootp server temporarily, powers on the 7100 CE using the initialization mode "Bootp & Store", waits for the Bootp cycle to be completed, closes the Bootp server and powers down the 7100 CE. Then, he selects the initialization mode "Using Stored" and powers on the 7100 CE again. From now on, he is able to establish the TCP/IP connection to the 7100 CE with the parameters obtained in that single Bootp cycle.

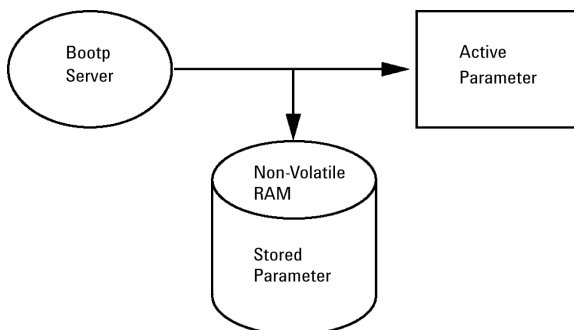
Figure 4 Bootp & Store (principle)

NOTE

Use the initialization mode "Bootp & Store" carefully, because writing to the non-volatile memory takes time. Therefore, if the 7100 CE is to obtain its parameters from a Bootp server every time it is powered on, the recommended initialization mode is "Bootp"!

Using Stored

When initialization mode "Using Stored" is selected, the parameters are taken from the non-volatile memory of the instrument. The TCP/IP connection will be established using these parameters. The parameters were configured previously by one of the described methods.

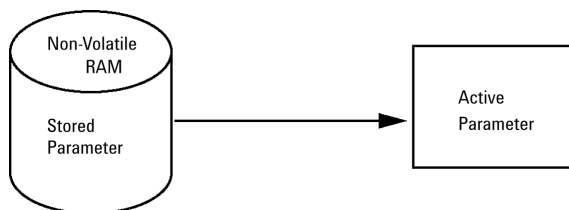
Figure 5 Using Stored (principle)

Using Default

When "Using Default" is selected, the factory default parameters are used instead. These parameters enable a TCP/IP connection to the LAN interface without further configuration, see [Table 9](#) on page 82.

A direct connection of a computer and the 7100 instrument is the entry point to change the stored parameters of the instrument. The default parameters of the 7100 instrument are permanently stored and cannot be changed. Using them, this allows you to establish a LAN connection at any time when the init mode "Using Default" is configured.

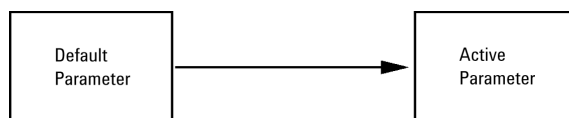
Figure 6 Using Default (principle)

NOTE

Using the default address in your local area network may result in network problems. Take care and change it to a valid address immediately unless you plan a direct LAN connection to stand-alone PC.

4 LAN Configuration

Initialization Mode Selection

Table 9 Using Default parameters

IP address:	192.168.254.11
Subnet mask:	255.255.255.0
Default gateway	Not specified

Since the default IP address is a so-called local address, it will not be routed by any network device. Thus, the PC and the CE instrument must be in the same subnet.

The user may open a Telnet session using the default IP address and change the parameters stored in the non-volatile memory of the CE instrument. He may then close the session, select the initialization mode "Using Stored", power-on again and establish the TCP/IP connection using the new parameters.

When the 7100 CE instrument is wired to the PC directly (for example using a cross-over cable or a local hub), that is separate from the local area network, the user may simply keep the default parameters to establish the TCP/IP connection.

NOTE

In the "Using Default" mode, the parameters stored in the memory of the 7100 Capillary Electrophoresis instrument are not cleared automatically. If not changed by the user, they are still available when you switch back to the mode "Using Stored".

Link Configuration Selection

The LAN interface supports 10 or 100 Mbps operation in full or half-duplex modes. In most cases, full-duplex is supported when the connecting network device - such as a network switch or hub - supports IEEE 802.3u auto-negotiation specifications.

When connecting to network devices that do not support auto-negotiation, the LAN interface will configure itself for 10 or 100-Mbps half-duplex operation.

For example, when connected to a non-negotiating 10-Mbps hub, the LAN interface will be automatically set to operate at 10-Mbps half-duplex.

If the CE instrument is not able to connect to the network through auto-negotiation, you can manually set the link operating mode using link configuration switches on the CE instrument.

Table 10 Link configuration switches

	SW3	SW4	SW5	Link configuration
	OFF	-	-	Speed and duplex mode determined by auto-negotiation
	ON	OFF	OFF	Manually set to 10 Mbps, half-duplex
	ON	OFF	ON	Manually set to 10 Mbps, full-duplex
	ON	ON	OFF	Manually set to 100 Mbps, half-duplex
	ON	ON	ON	Manually set to 100 Mbps, full-duplex

Storing the Settings Permanently with Bootp

If you want to change parameters of the CE instrument using Bootp follow the instructions below.

- 1 Turn off the CE instrument.
- 2 Change the CE instrument's settings for the configuration switch to "*Bootp & Store*" mode, see [Table 8](#) on page 79.
- 3 Start the Agilent Bootp Service (see "[Agilent Bootp Service](#)" on page 89) and open its window.
- 4 If required, modify the parameters for the CE instrument according to your needs using the existing configuration.
- 5 Press OK to exit the Bootp manager.
- 6 Now turn on the CE instrument and view the Bootp server window. After some time the Agilent Bootp Service will display the request from the LAN interface. The parameters are now stored permanently in the non-volatile memory of the CE instrument.
- 7 Close the Agilent Bootp Service and turn off the CE instrument.
- 8 Change the settings of the CE instrument's configuration switch to "*Using Stored*" mode, see [Table 8](#) on page 79.
- 9 Power cycle the CE instrument. The CE instrument can be accessed now via LAN without the Agilent Bootp Service, refer to "[Agilent Bootp Service](#)" on page 89.

Manual Configuration

Manual configuration only alters the set of parameters stored in the non-volatile memory of the CE instrument. It never affects the currently active parameters. Therefore, manual configuration can be done at any time. A power cycle is mandatory to make the stored parameters become the active parameters, if the initialization mode selection switches permit this.

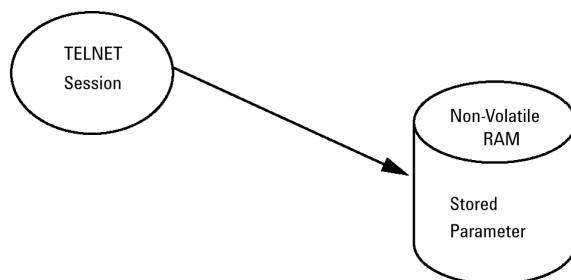
Figure 7 Manual configuration (principle)

With Telnet

Whenever a TCP/IP connection to the 7100 CE instrument is possible (TCP/IP parameters set by any method), the parameters may be altered by opening a Telnet session. Please refer to “[Using Default](#)” on page 81 for establishing an initial LAN connection.

- 1 Open the system (DOS) prompt window by clicking the Windows **START** button and select "**Run...**". Type "cmd" and press OK.
- 2 Type the following after the system (DOS) prompt:

- c:\>telnet <IP address>
- or
- c:\>telnet <host name>

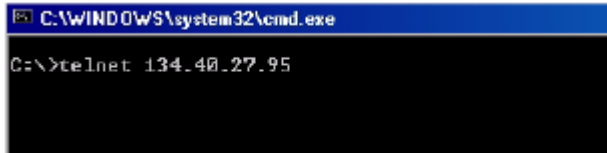
Figure 8 Telnet - starting a session

where <IP address> may contain the assigned address from a Bootp cycle, or the default IP address (see “[Configuration Switch](#)” on page 78).

If the connection is established successfully, the instrument responds with the following:

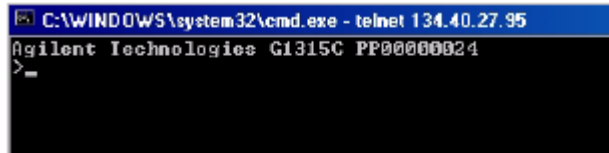
Figure 9 A connection to the instrument has been made

3 Type

?

and press **[Enter]** to see the available commands.

```

C:\WINDOWS\system32\cmd.exe - telnet 134.40.27.95
Agilent Technologies G1315C PF00000024
>?
command syntax          description
-----
?                        display help info
/                        display current LAN settings
ip <x.x.x.x>             set IP Address
sm <x.x.x.x>            set Subnet Mask
gw <x.x.x.x>            set Default Gateway
exit                    exit shell
>
  
```

Figure 10 Telnet commands

Table 11 Telnet commands

Value	Description
?	Displays syntax and descriptions of commands
/	Displays current LAN settings
ip <x.x.x.x>	Sets new IP address
sm <x.x.x.x>	Sets new subnet mask
gw <x.x.x.x>	Sets new default gateway
exit	Exits shell and saves all changes

4 To change a parameter follows the style:

- parameter value, for example:
ip 134.40.27.230

Then press **[Enter]**; parameter refers to the configuration parameter you are defining, and value refers to the definitions you are assigning to that parameter. Each parameter entry is followed by a carriage return.

4 LAN Configuration

Manual Configuration

- 5 Use “/” and press [Enter] to list the current settings.

```
C:\WINDOWS\system32\cmd.exe - telnet 134.40.27.95
>/
LAN Status Page
-----
MAC Address   : 0030D30A0038
Init Mode    : Using Stored
-----
TCP/IP Properties
- active -
IP Address   : 134.40.27.95
Subnet Mask  : 255.255.248.0
Def. Gateway : 134.40.24.1
-----
TCP/IP Status : Ready
-----
Controllers  : no connections
>=
```

Information about the LAN interface
MAC address, initialization mode
Initialization mode is Using Stored
active TCP/IP settings

TCP/IP status - here ready
connected to a PC with controller software
(for example Agilent ChemStation), here not
connected

Figure 11 Telnet - current settings in "Using Stored" mode

- 6 Change the IP address (in this example 134.40.27.99) and type “/” to list the current settings.

NOTE

If the initialization mode switch is changed now to “Using Stored” mode, the instrument will take over the stored settings when the instrument is re-booted. In the example above, it would be 134.40.27.99.

Agilent Bootp Service

Agilent Bootp Service provides central administration of IP addresses for Agilent CE instruments in a LAN. The service runs on the instrument LAN PC, which must be running a TCP/IP network protocol and cannot run a DHCP server.

When an instrument is powered on, a LAN interface in the instrument sends a request for an IP address or host name and provides its hardware address as an identifier. The request may continue for up to 5 minutes. The Agilent Bootp Service answers this request and passes a previously defined IP address and host name associated with the hardware address to the requesting instrument.


When the instrument receives its IP address and host name, it stops sending the request. It maintains the IP address as long as it is powered on. Powering down the instrument causes it to lose its IP address, so the Agilent Bootp Service must be started. Since the Agilent Bootp Service runs in the background, the instrument will receive its IP address once it is powered up.

Detailed documentation on installation and the usage of the Agilent Bootp Service can be found on your ChemStation installation DVD under **\Support\Manuals\Installation**.

Location of the Media Access Control (MAC) Address

The MAC or hardware address of the LAN interfaces is a world-wide, unique identifier. No other network device will have the same hardware address. The MAC address is physically located on the electronic main board in the rear box of the instrument. It is not accessible by the user.

The MAC can be reviewed from a connected instrument by accessing it via a Telnet session as described earlier. Please follow the instructions on Manual Configuration, Using Telnet (“[With Telnet](#)” on page 86) and see [Figure 11](#) on page 88 for the location of the MAC address as the second line in the LAN status page.

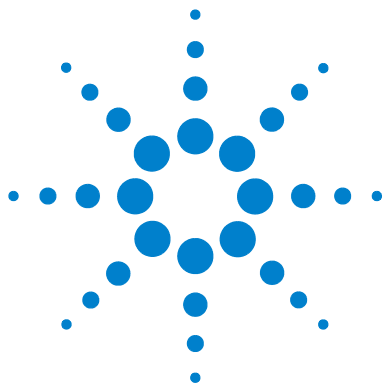5 Using the Agilent 7100 Capillary Electrophoresis System

System Definition	92
Controlling the Agilent CE Instrument	93
Working with the CE Diagram	98
How to Prepare your Agilent CE Instrument for Analysis	105
Capillary Electrophoresis, CE	106
Capillary Electrophoresis, Plus High Pressure CE+p	157
Capillary Electrochromatography, CEC	169

This chapter provides information on how to set up the CE instrument for an analysis and explains the basic settings.

System Definition

After your Agilent 7100 Capillary Electrophoresis System has been properly installed (see “[Installing the Instrument](#)” on page 67) you can begin exploring the system. Your Agilent 7100 Capillary Electrophoresis System comprises the CE instrument, the ChemStation and the Lab Advisor/Instrument Utilities software. The CE instrument doesn't have a keyboard itself. It is completely controlled by the ChemStation through the CE and DAD diagram.

For detailed description of data evaluation see the manuals “Understanding Your ChemStation” and “Agilent ChemStation for CE System, Installing and Understanding”

Controlling the Agilent CE Instrument

Before You Start the Agilent CE Instrument

Use the following as a checklist before you start the Agilent CE instrument.

Make sure that:

- the replenishment bottles are installed and properly tightened,
- the cassette (with a capillary installed) is inserted in the system,
- the cassette door and the tray door are closed, and
- you do not need to put the vials into the tray before you start the Agilent CE instrument

Starting Up the Agilent CE Instrument

The power switch of the instrument is located on the left front corner of the instrument.

Press this switch to turn on the line power.

After the instrument is started, the fans and the pump start working. The instrument starts to build-up internal pressure (approximately 950 mbar) first, followed by a vacuum. The pump becomes idle and switches on only occasionally once the pressure and vacuum setpoints are reached.

During initialization the instrument tests different functions and adjustments. The tray as well as all three lift stations are moved.

The UV lamp will switch on during instrument startup based on parameters that can be defined in the DAD control section, see “[Diode Array Detector](#)” on page 199. Otherwise, the lamp can be switched on from the DAD diagram of the ChemStation.

The meaning of the LED colors on the front cover are explained in the following table.

Table 12 Status LED color

Color	Meaning
No color	Instrument is ready
Green	Analysis running, do not use instrument with the exception of manual operation of the tray. Some fields in DAD and CE diagram are shaded gray.
Yellow	Not-ready condition. The run cannot be started yet. When you start, the LED will come on after a certain time. Click the Agilent CE System status bar in the ChemStation to see what is not ready (for example temperature).
Red	Error condition. There was or is currently a problem with the instrument or communication to the PC, which requires your action or an acknowledgement. To get more information, connect to the instrument with the Lab Advisor or ChemStation software. Check the logbook.

NOTE

A blinking red LED indicates the instrument is in a resident mode. This means only a core of the firmware is active and the instrument has limited functionality. The Lab Advisor software will be able to connect to the instrument under these circumstances. An instrument restart by switching the power off and on again boots with complete firmware loaded. During a main firmware download (see “[Firmware](#)” on page 196), the instrument shows a red blinking LED.

Starting Your Computer and the Agilent ChemStation Software

After you have switched on your computer and it has booted up, start the Agilent ChemStation by double clicking the **Instrument Online** icon in the Agilent ChemStation menu.

NOTE

In order to control the 7100 Capillary Electrophoresis instrument, the ChemStation revision must be C.01.01, MassHunter Acquisition for TOF/QTOF B.05.01, MassHunter Acquisition for QQQ B.06.00 or higher.

NOTE

We recommend you to install the latest Lab Advisor software revision (B.02.09 or higher).

If you have problems starting the Agilent ChemStation, check that the software has been properly installed. During startup the ChemStation might ask you to update the connection settings, see [Figure 12](#). You might suppress this message in the future. If you want to update the instrument configuration, select **Yes**. Typically, an instrument configuration must be updated for first time use of the software, changes in the IP address or changes in the mode of usage.

NOTE

Please refer to the ChemStation documentation and online help for details on installation and configuration of the 7100 CE instrument.

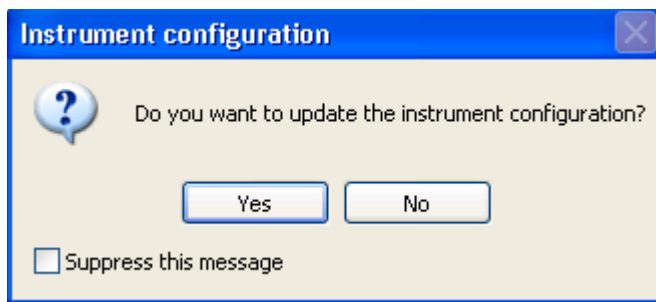
Figure 12 Instrument configuration during ChemStation startup

Starting the Agilent ChemStation Software the First Time

Starting the Agilent ChemStation software the first time will lead to the following message:

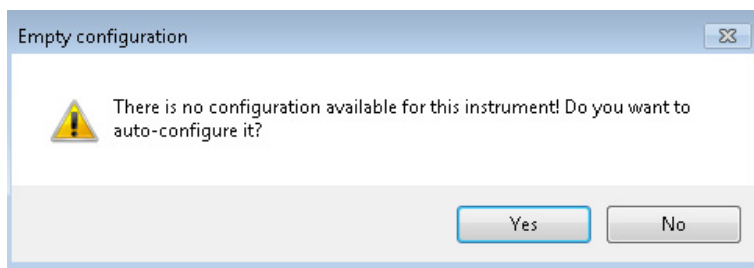
Figure 13 Configuration at first start of ChemStation

Please select **Yes**. The following dialog will allow you to configure the instrument. Please provide an IP address or host name in the subsequent dialog. Check the respective radio button and type in either the IP address or the host name of your instrument.

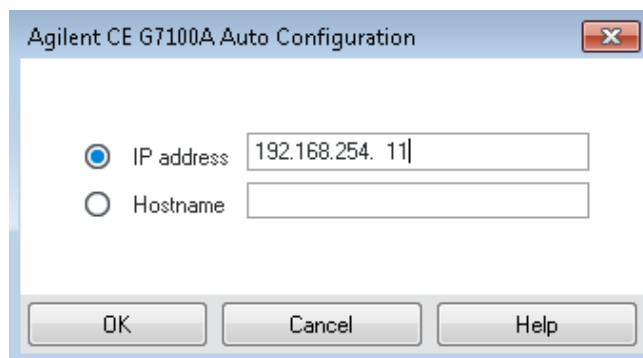
Figure 14 Request of IP address or host name

The auto-configuration process will detect an instrument at the IP address or host name you provide. You can confirm the instruments identity with the serial number.

Please select "Configure..." if you would like to connect to a different instrument.

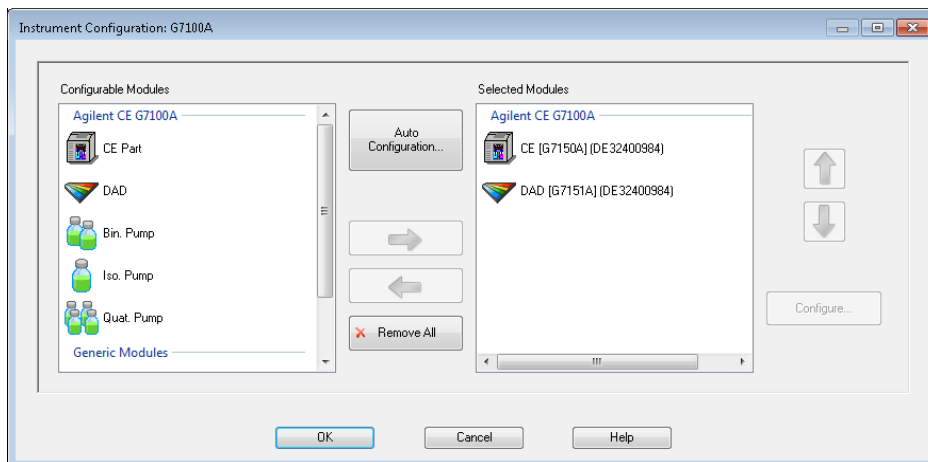
Figure 15 Configuration menu

NOTE

Please refer to “[LAN Configuration](#)” on page 75 for the default IP address or manual configuration of an individual IP address.

After you clicked the Launch in the OpenLAB control panel, the software loads the last used method. The software boots in the exact last status.

Do not use the instrument or the graphical user interface during initialization. Wait until the system has reached the ready state.

If there is a vial in tray position 50 (reserved position), the system asks you to remove that vial before you proceed.

Duplicate vials are put in the tray and the respective position is brought to the front. Duplicate vials may occur if vials from the tray are loaded in a lifter and later on another vial is loaded in the same tray position.

If there are no bottles (electrolyte or waste) in the replenishment system or they are not properly tightened, the error LED (red) will come on because no pressure can be built up. Check the logbook for more information.

Working with the CE Diagram

When you start the Agilent ChemStation, the CE and DAD diagram appears by default as the main screen. The CE diagram displays the Agilent CE instrument showing the current status of the system as a whole. Each icon represents one component in the system.

Common Software Toolbar

The software tool bars in the Agilent ChemStation offer a group of commonly used functions. The software tools bars visible in the Method and Run Control View are shown in [Figure 17](#) on page 100. To change values for a component, you must choose the appropriate icon.

This will bring up the menu. The menu may comprise sets of items; for changing parameters in your method, for control of the instrument, and for help.

The tool bars associated with the Data Analysis View and Report Layout view look similar and are described in the online help and the ChemStation documentation.

For more assistance, click **Help** in the menu bar.

Software Toolbar

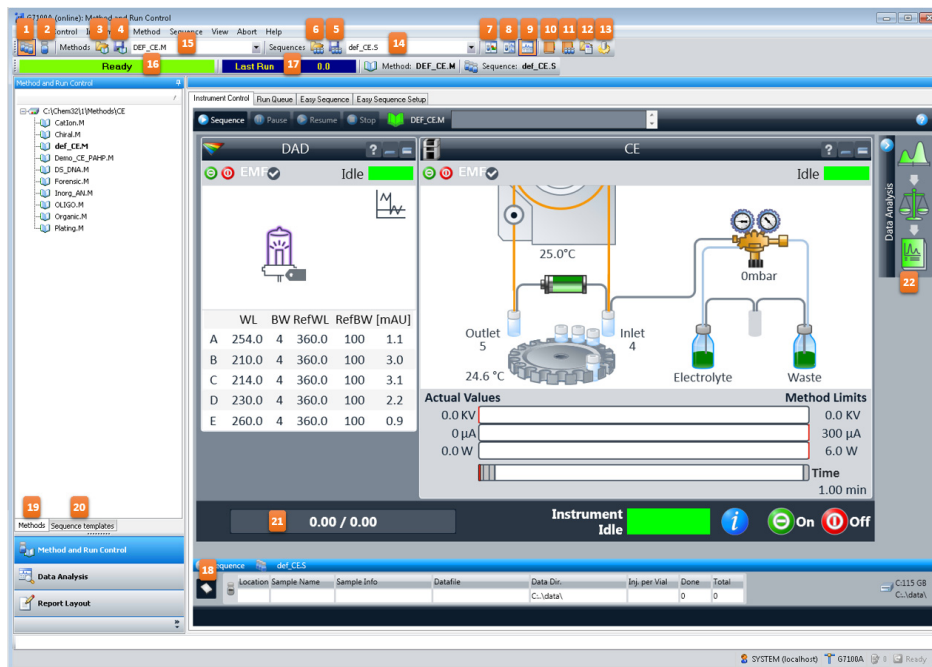
Figure 16 Components of the common software toolbar of the ChemStation.

- | | | |
|--|--|---|
| 1 Work with sequences | 9 Change the layout to show the system diagram | 17 Display of elapsed and overall runtime |
| 2 Work with single vial analysis | 10 Display the current logbook | 18 Display sample information for sequences or single vials |
| 3 Load method | 11 Display the current sequence logbook | 19 Access method templates |
| 4 Store method | 12 View latest report file(s) | 20 Access sequence templates |
| 5 Store sequence | 13 Open the Preferences dialog | 21 Display of elapsed and overall runtime |
| 6 Load sequence | 14 Name of current sequence and quick access to recent sequences | 22 Access integration, calibration and reporting settings |
| 7 Change the layout to show the sampling area and system diagram | 15 Name of current method and quick access to recent methods | |
| 8 Change the layout to show two online plots | 16 Status bar | |

Graphical User Interface Explanation

Components of the Agilent CE instrument in the CE diagram (synonym dashboard) are shown in [Figure 17](#). Each component in the instrument is represented by an icon in the CE diagram or DAD diagram.

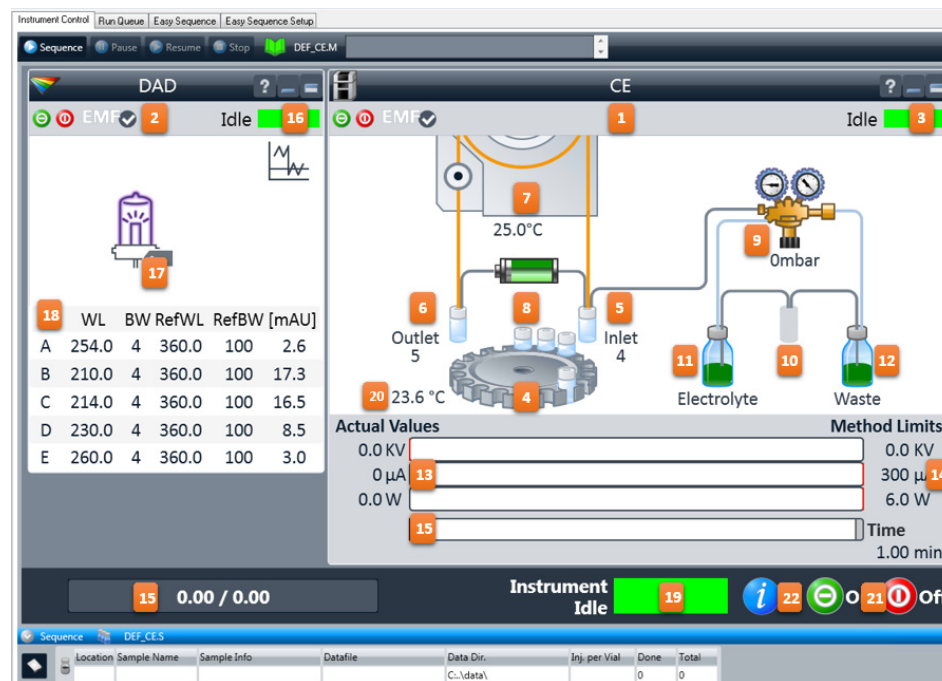
Figure 17 The CE and DAD diagram represents each component in the instrument

- | | | |
|-----------------------------------|--|-------------------------------------|
| 1 CE control diagram | 9 Internal pressure | 17 Lamp & Lamp TAG information |
| 2 DAD control diagram | 10 Replenishment station | 18 Detector wavelengths/spectra |
| 3 CE status | 11 Electrolyte bottle / pressure storage | 19 Instrument status |
| 4 Tray | 12 Waste bottle / vacuum storage | 20 Tray temperature |
| 5 Inlet lift / loaded vial | 13 Current electrical values | 21 Device control (ready/off) |
| 6 Outlet lift / loaded vial | 14 Electrical parameter from method | 22 Display instrument configuration |
| 7 Cassette temperature | 15 Acquisition | |
| 8 High voltage power supply state | 16 DAD status | |

The current analysis mode is displayed in the top line of the CE diagram. The analysis modes that can be selected under instrument configuration are CE, CE+p, CEC or CE-MS.

For configuration without a DAD detector, for example with external detectors, the DAD diagram will not be displayed. For configuration with an additional pump, for example CE-MS, an additional pump diagram will be displayed.

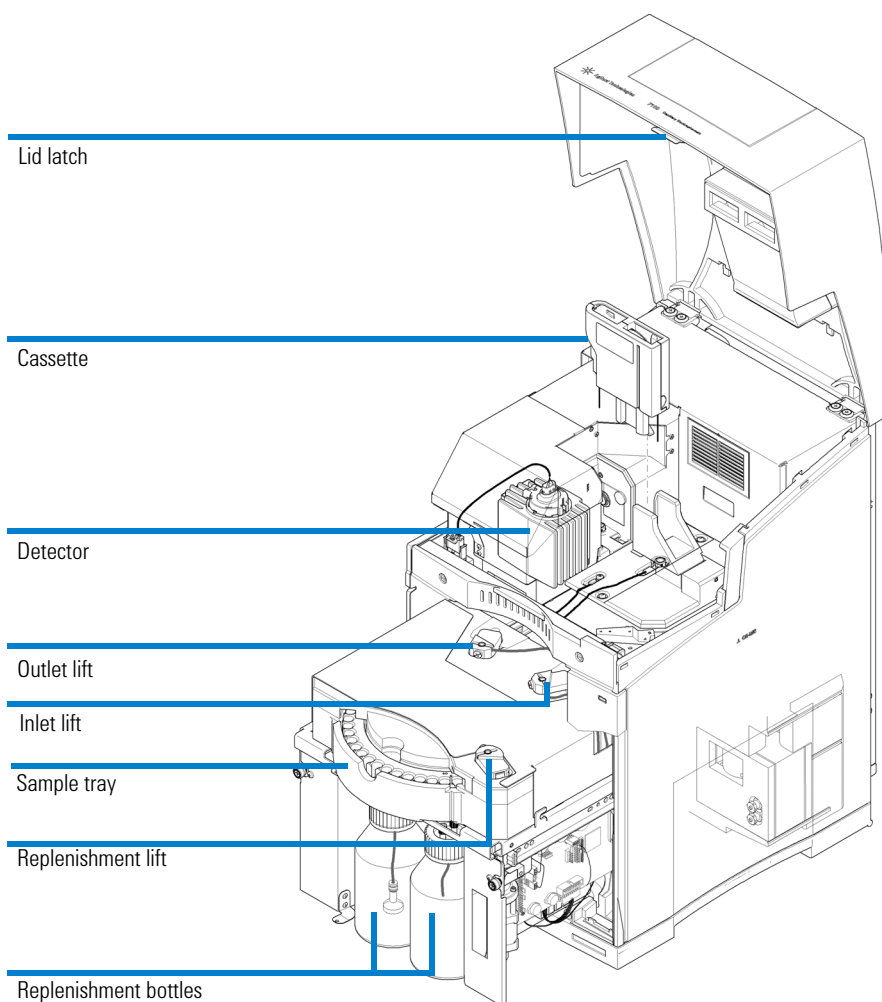
Figure 18 Corresponding hardware components

Menus in the CE and DAD Diagram

The CE and DAD diagram offers access to directly control several instrument components, as well as access to method parameters. Right-clicking in the background of each diagram provides access to the context menus (Figure 19).

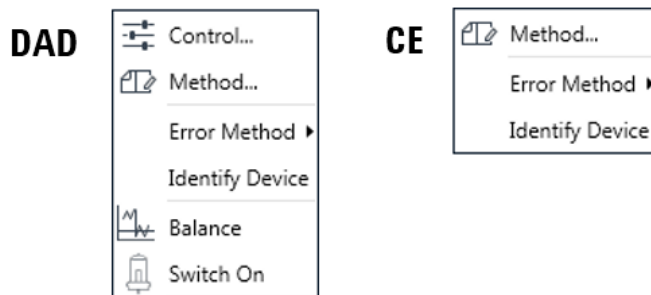
Figure 19 Menus in the DAD and CE diagram

To trigger actions for an instrument component, select the appropriate icon in the diagram. A context menu including possible actions, such as switching on the lamp, appears. Any instrument action triggered through such a menu does not change the current method.

To change values for a component, select the appropriate icon. A menu with three general items and parameter specific for the component appears.

Method

A click here requests the current method from the instrument and opens the CE **Method Setup** dialog box for the CE part of the current method or the respective DAD part. To set up a complete method, refer to “[How to Use Your Agilent CE Instrument for Capillary Electrophoresis Separations](#)” on page 106. If closed with **OK** the (possibly) modified method is downloaded to the device. If changes have been made an asterisk is added to the method status. The method status is represented by a book icon in the top line of the dashboard.

Direct Control of Instrument Components

The actions are performed immediately and the changes are reflected in the CE diagram. Such actions include, for example:

- Switching the lamp on.
- Controlling the sample tray.
- Lowering the lift stations to remove the capillary cassette.
- Releasing the pressure to change the replenishment bottles.

Working with these items does not change the method currently loaded.

Set Error Method

Set Error Method enables you to store the current parameters. These parameters are applied to the instrument when an error occurs. It ensures that the instrument shuts down in a controlled manner if ChemStation control is discontinued for any reason.

Identify Device

The command **Identify Device** causes the LED on the front cover to blink alternating in green and yellow for 4 seconds.

System Vialtable

The vialtable is a graphical tool which facilitates the management of vials in the tray. It can be used as an intelligent notepad, which exchanges its information with methods and sequences.

Some features are:

- If a vial is defined in the vialtable this definition (name of vial) will be used whenever the vial is referenced (method editing, method print).
- Consistency checks can be done in order to verify if there are vial conflicts between the method and sequence usage of vials.
- Documentation of vial positions and contents.

How to Prepare your Agilent CE Instrument for Analysis

To prepare the Agilent CE instrument for the analysis you need to do the following tasks:

- Start the Agilent CE instrument.
- Start the computer.
- Start the Agilent ChemStation software/CE user interface.
- Allow the lamp to warm up (approximately 1 hour).
- Prepare vials for sample and buffer.
- Insert a capillary into the alignment interface.
- Insert the capillary into the capillary cassette.
- Insert the capillary cassette into the system.
- Prepare the replenishment system (optional).

After you have completed all these tasks, you can start an Agilent CE method.

A CE method can be created in an online (CE instrument is on and connected) or offline installation of the Agilent ChemStation.

The detector lamp can be switched on automatically and unattended by scheduling this from the detector control menu. Please see the options for the detector under “[Diode Array Detector](#)” on page 199.

The detector lamp can be switched on automatically with each instrument start. Select the respective setting from the detector control menu (see “[Diode Array Detector](#)” on page 199).

Capillary Electrophoresis, CE

How to Use Your Agilent CE Instrument for Capillary Electrophoresis Separations

This chapter introduces method parameters for applications running in CE mode. This mode is the most suitable for applications in the field of capillary zone electrophoresis (CZE) and micellar electrokinetic chromatography (MEKC). The current CE mode for your instrument can be selected under **Instrument > more CE > Instrument configuration...** . The current mode is visible in the second line of the CE control diagram. For detailed information, see [“Working with the CE Diagram”](#) on page 98 and [“Graphical User Interface Explanation”](#) on page 100.

Agilent CE Method Parameters

This section describes the parameters of an Agilent CE method for data acquisition and their meanings. Example sets of parameters for a method to analyze the test sample are in [“Setting Up a Test Sample Run”](#) on page 272. After you have completely set up the Agilent CE method and supplied the sample information, you can run your method. Choose **Run Method** from the Run control menu.

Overview

A convenient way to access the parameters of an Agilent CE method is the item **Setup Instrument Method** in the **Instrument** menu. This displays the **CE Method Setup** dialog box. Tabs with parameters for the DAD and CE part of a method can be edited on this screen. The left-hand side shows the most important parameters. The right-hand side offers several pages that can be expanded subsequently for editing.

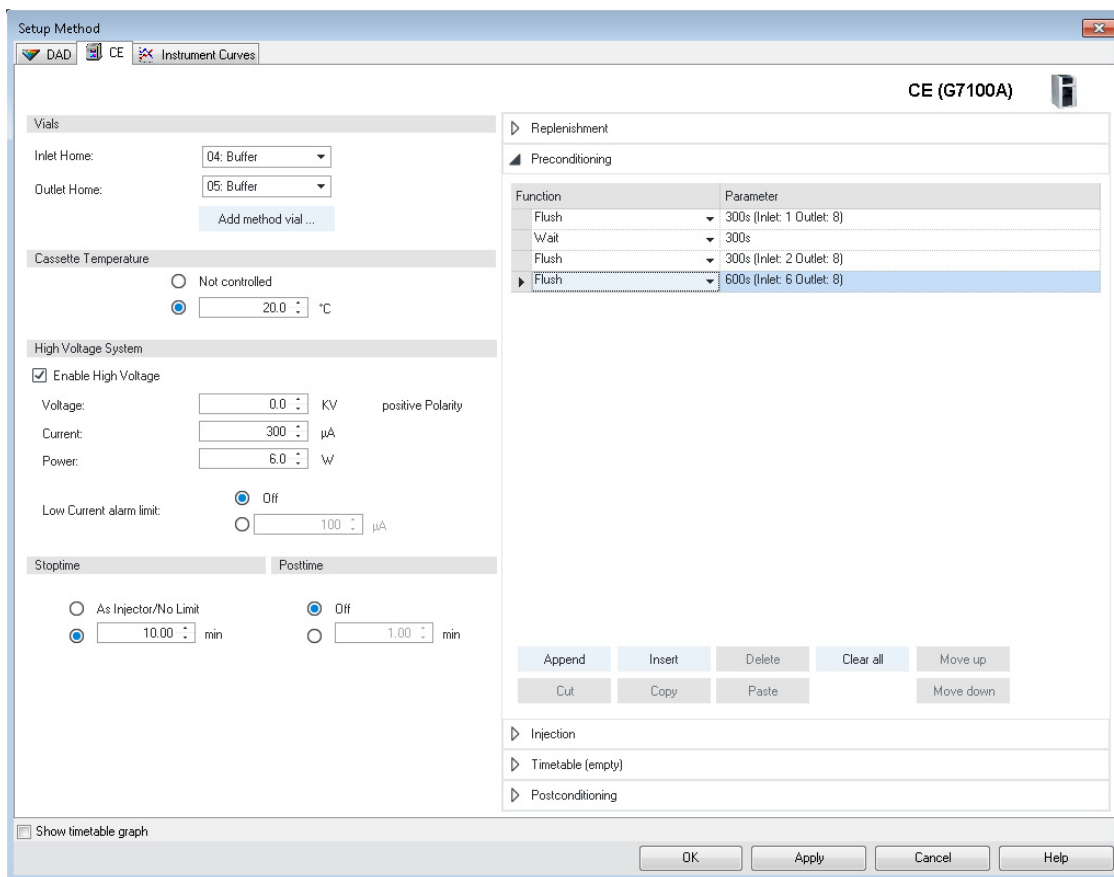
Figure 20 CE Method Setup

5 Using the Agilent 7100 Capillary Electrophoresis System Capillary Electrophoresis, CE

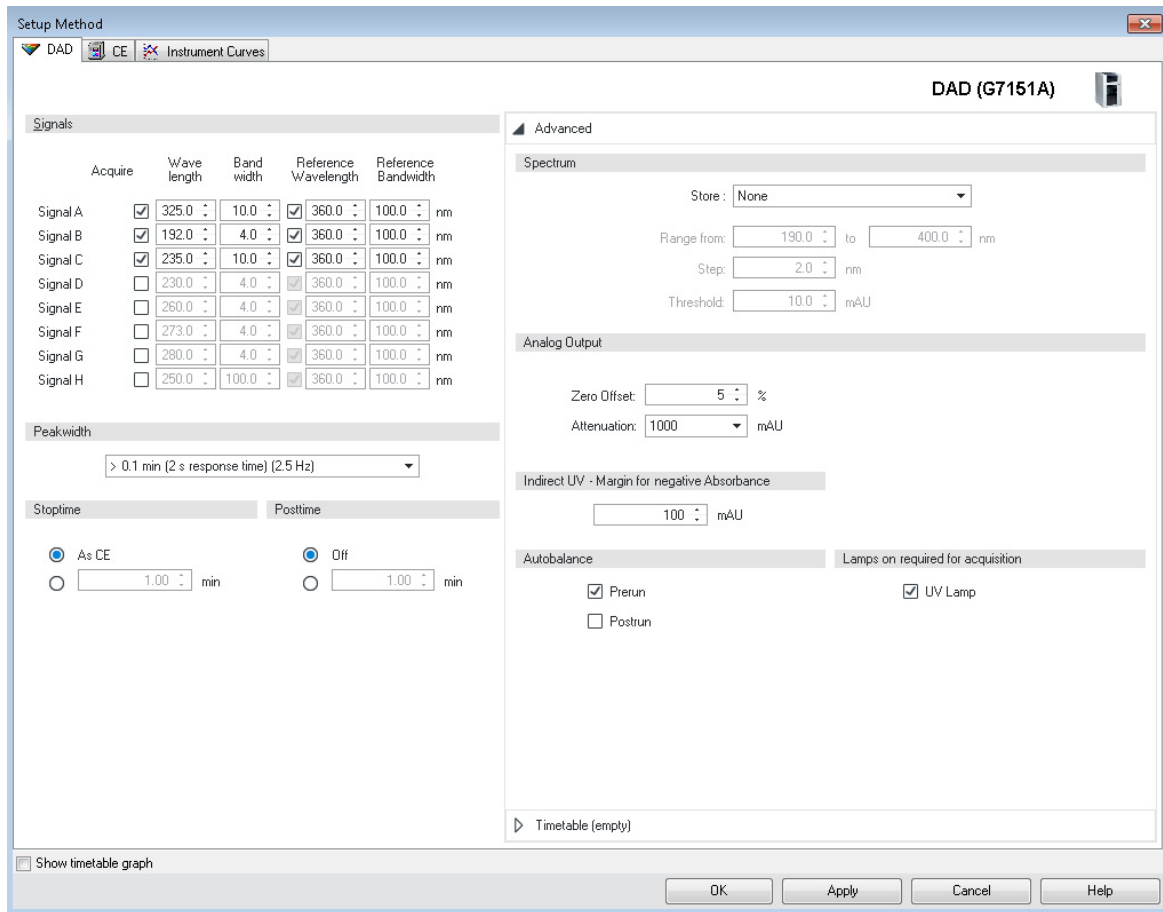
Figure 21 DAD Method Setup

Selecting **OK** transfers the parameters to the instruments and closes the dialog. Choosing **Apply** only transfers the current parameter to the instrument. Save the edited method, if appropriate.

If you do not save the method to the hard disk, this changed method is executed when choosing **Run Method**. When you refer to a method in a sequence, the method is always loaded from the disk. So if you did not save your changes using **Save Method**, the original method from the hard disk is used.

Method parameters are also available if **Method/Edit Entire Method...** is chosen. This dialog guides you through further parameters associated with a method such as

- Method Information
- Instrument / Acquisition
- Data Analysis
- Run Time Checklist

The book icon in the software toolbar is a further alternate entry point for method editing. When you click the icon, you can access menu items to set method parameters.

Vials

You can program a method to use certain vials as the inlet or outlet home vial. During electrophoresis, the electrodes and the capillary are in these vials. Vial numbers 1 to 48 or user vials (User 1 to User 10) can be selected. The selection can be done from the pull-down menu or by typing in the number with the keyboard. Clicking **Add Method Vial** allows you to enter vial names which will be visible in the vialtable.

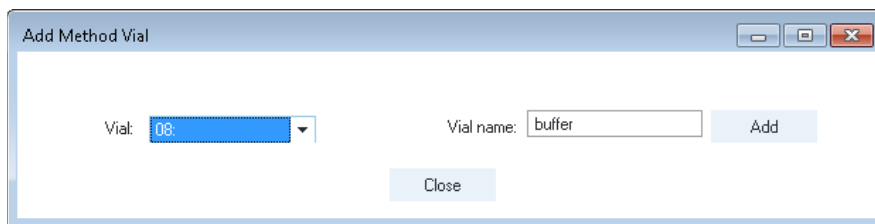
Figure 22 Add Method Vial

User vials are numbered from 1 to 10. User vials are placeholders in the method which are resolved by the sequence table at a later point in time. Usage of user vials is helpful for method optimization. Various buffer compositions can be tested with one common method. Creation of methods with variations of Inlet or outlet home vials can be omitted. Usage of user vials is also helpful to change buffers in long sequences to avoid over utilization of the buffer.

NOTE

Vial numbers below 10 must be typed in with a preceding 0, for example 02 for a vial in tray location number 2, since the vials are numbered with two digits.

Cassette Temperature

The cassette temperature can either be set to not be controlled or to a certain setpoint. This setpoint must be within the instrument's specified range (see "Physical Specifications" on page 21).

The lowest applicable temperature is 10 °C below ambient temperature. The highest temperature is 60 °C.

The instrument is ready after the cassette temperature reaches the setpoint and stabilizes. A run can be started.

The home values are activated at five points in your analysis:

- When a method is loaded.
- When you activate **Run Method** or **Run Sequence** from the **RunControl** menu.
- After injection is completed.
- After the stoptime has elapsed.
- After the entire analysis (including postconditioning) is completed.

NOTE

Keep the cassette temperature above the dew point or decrease the ambient humidity.

The instrument is specified to be operated at non-condensing conditions which is at humidity below 80% at 31 °C (87.8 °F). Please be aware: if the temperature of the cassette is below the dew point condensation might take place. Subsequently water condensate from the ambient might lead to arcing and damage of the cassette or instrument.

High Voltage System

The following parameters for the power supply can be controlled using the CE **Method Setup** screen in the Agilent ChemStation software:

- Enable High Voltage
- Voltage
- Current
- Power

In addition you can set a Lower Alarm Limit for Current.

NOTE

In order to directly apply HV from the dialog associated with the HV icon (a battery) in the CE dashboard, the active instrument method must have **Enable High Voltage** check marked.

NOTE

The maximum stored energy in the CE power-supply is below 200 mJ at 30 kV. Due to this and the fixed current limit of 300 μ A the high voltage applied is not defined as hazardous voltage.

Positive voltage settings will configure the electrode at the inlet vial with positive values (positive polarity). Positive polarity is regarded as the standard setting.

You can reverse the polarity to negative by entering negative values for high voltage. This means that a negative voltage is applied at the inlet electrode.

The outlet electrode is always grounded to ensure that the potential at the point of detection is close to ground. Even with negative polarity, the inlet electrode is the point where the voltage is applied and the outlet electrode is close to ground.

Although the outlet electrode is always at the ground potential, you should make sure that the electrode is not short-circuited to ground. The electrode is connected to the power supply to measure the current flowing through the capillary. If the ground electrode is short-circuited to ground, all current flowing through the capillary is signaled as leakage current.

NOTE

It is good CE practice to ramp up voltage instead of switching voltage suddenly to high values. This is regarded as gentle treatment for samples which might otherwise be affected by the field strength, such as proteins. A linear voltage ramp from 0 to 30 kV can be programed by starting with 0 kV and setting the voltage in the timetable to for example 0.2 minute.

Voltage, Current and Power

Voltage, current and power are related to each other by the resistance of your capillary/buffer system. You can set values for each of these three parameters. The individual values are treated as limits. The limit that is reached first applies. You can only enter absolute values for current and power.

If you only want to control the voltage and not the current and power, you can set those two to the system limit (typically: 300 μ A, 6 W) and the system will permit the maximum value.

Lower Alarm Limit for Current

When using constant voltage mode, the current is normally the indicator of the stability of the system. A rapid current decrease or drop can indicate instability in the system. By setting a lower alarm limit you can define when you want to be warned of such instable conditions during the run.

Instable conditions can be caused by:

- No buffer in the capillary
- Air bubble in the capillary
- Clogged capillary
- Broken capillary

Current values are dependent on the buffer concentration, the type of buffer used, the inner diameter and length of the capillary, and the temperature. When running the Example Method for the Test Sample (see [“Setting Up a Test Sample Run”](#) on page 272) normal current values are below 50 μ A.

If you set the lower alarm limit and the it is not exceeded 30 seconds after start of the run, the run is stopped. When the lower alarm limit is reached during the run part of the analysis (when the HV is applied), your current run is stopped. During a sequence the system continues with the next analysis.

Do not use the lower alarm limit or program the time if changing the inlet or outlet vial during the run (for example for fraction collection). Otherwise, the system will stop when changing the vial because the lower alarm limit is reached.

Posttime and Stoptime

The stoptime defines the time for your separation, also referred to as the run part of the analysis. It does not include replenishment, preconditioning, injection and postconditioning. After the stoptime has elapsed, the system changes from the run state to postconditioning or postrun (if defined) and then to the ready state.

The posttime is normally used for equilibration. It defines the time after the analysis, that is after the stoptime has elapsed and data analysis is completed. During the posttime the instrument is in the postrun state. After both stoptime and posttime have elapsed the instrument completes postconditioning and is then ready for the next analysis.

Replenishment

During multiple analyses, the composition of the buffer can change, for example the pH due to electrolysis. A simple way to refresh the buffer is to access different buffer vials in the tray which might be “user vials” in the method that are assigned to a certain tray position in the sequence table. This is more useful, for example, during method development.

Another way to change the buffer is replenishment. Replenishment can be programmed via the topmost table in the right side of the CE **Method Setup** screen. It is done either in parallel or serial to the subsequent conditioning operation. The associated timetable can be used to empty, fill or replenish vials to a certain height.

The replenishment system automatically empties and refills the buffer vials in the tray with new buffer from the electrolyte bottle. Therefore, the number of tray positions occupied by home vials is reduced and a maximum of sample vials can be loaded. Typically replenishment is done after every 3 or 4 runs. Depending on the stability of the running buffer, it is sometimes necessary to do it before each run. If you want to use the replenishment system, prepare it as described in [“Preparing the Replenishment System if Needed”](#) on page 125.

NOTE

Using parallel replenishment, you can decrease the total time needed for your analysis, as replenishment and preconditioning are done at the same time. When using this feature, you have to make sure that vials used in conditioning are not replenished at the same time, otherwise you will have a vial conflict.

NOTE

The time for a replenishment step can vary from vial to vial.

When using buffers containing detergent (for example SDS), replenishment may not fulfill your needs. The limit for SDS concentration is 70 mM. Do not exceed this concentration. Add some organic liquid to the waste bottle so that bubbles do not form. Also see [“Problems with the Replenishment System and Buffers Containing SDS”](#) on page 188.

Preconditioning

In order to have stable conditions for your analysis, it is important to condition capillaries.

Depending on the separation mode used it can comprise:

- conditioning the capillary with conditioning agents, and
- equilibrating the capillary with buffer.

These conditioning steps can be programmed using an expandable table in the right side of the CE **Method Setup** screen. The associated timetable can be used to apply pressure, apply voltage or flush between vials. Inlet and outlet vials can be loaded and wait time can be programmed here. If suitable, such steps can also be done as a postconditioning operation ([“Preconditioning and Postconditioning”](#) on page 164).

NOTE

Condition capillaries according to their specifications. Fused silica capillaries are typically rinsed with alkaline or acidic solutions for conditioning. Please be aware that most coated capillaries will be irreversibly damaged by this.

NOTE

To minimize contamination of buffer vials with conditioning solutions (such as NaOH) which might stick to the outside of the capillary or electrode, it is recommended to dip the electrode in an additional step in a large dedicated vial filled with buffer prior to letting the instrument load the actual inlet and outlet home vials with run buffer. This precaution can be programmed into the timetable by a combination of the commands “load vial” and “wait”.

Wash Inlet Electrode (New Command to Avoid Carry-over in Critical Applications)

During precondition and postconditioning a command for emptying the space between capillary and electrode can be executed. The command is dedicated for applications with aqueous buffers and consists of application of slight vacuum to the inlet vial while lowering the inlet vial. This leads to opening of an internal sealing of the electrode and an airstream inside the wash vial at the inlet lift. Liquid stored in the space between capillary and electrode is removed. It is recommended to perform this prior to a flush step with run buffer in case of carry-over critical applications. The wash inlet electrode command is not available for the outlet electrode. At the inlet lift a dedicated wash vial should be present. Using the Wash Inlet Electrode commands leads to injection of a small volume on the outlet side (equaling approximately 50 mbar x 3 s for a 50 µm capillary). Therefore, position at the outlet side a vial with run buffer or any other suitable run buffer vial. Empty vials at the outlet lead to air in the capillaries which needs to be flushed out again.

NOTE

Usage of short electrodes (part number **G7100-60033**) in combination with pressure injections also reduces this type of take over greatly (insert #reference to injection section). Short electrodes are shorter by 5 mm, sample is filled only at low level and so the electrode does not touch the sample.

Injection

You have several choices for CE injection:

- Hydrodynamic injection (by pressure or vacuum)
- Electrokinetic injection (applying a voltage, current or power)
- Using an injection program

The injection parameters can be programmed using an expandable table in the right side of the CE **Method Setup** screen. The associated timetable can be used to apply pressure, apply voltage or flush between vials. Inlet and outlet vials can be loaded and wait time can be programmed here. After the injection has finished, the analysis starts with recording detector signals to result file.

Hydrodynamic Injection

With hydrodynamic injection, the inlet buffer reservoir is replaced with the sample vial. A pressure is applied for a certain time to introduce the sample in the capillary. The system constantly controls the pressure and corrects for the rise time effects of valves and the compressibility of air. When pressure is applied, the pressure to the sample vial is increased gradually to its setpoint after which the pressure decreases gradually to approximately a fifth. Then a correction time is inserted after which the pressure decreases gradually to atmospheric pressure. This results in accurate and reproducible injection, as well as exceptional injection linearity.

Injection by pressure is the most frequently used injection technique. There are no differences in injection concentration for molecules with different mobilities as in electrokinetic injection.

With viscous buffer system the flush command can be used. However, injection of sample to the capillary at this high pressure is not monitored by the self-correcting injection algorithm. This sample application to the capillary is therefore less reproducible. To compensate for this, internal standards might be useful in quantitative analysis.

Electrokinetic Injection

For electrokinetic injection the inlet buffer vial is replaced with the sample vial. A voltage, current or power is applied for a certain time that causes the sample to migrate into the capillary.

This injection technique is often used for capillaries filled with fixed or cross-linked gels or other high-viscosity material, where injection by pressure is not applicable.

During electrokinetic injection, molecules with different mobilities will be depleted differently from the sample solution, therefore gradual changes in analyte composition might occur as a consequence of multiple injections.

Different salt contents in the sample matrix will influence the injection efficiency of the analytes. Variation in the matrix composition might therefore bias the quantitative analysis and, in the worst case, hamper analyte introduction.

Difference Between Hydrodynamic and Electrokinetic Injection

To do electrokinetic injection, the electrode must touch the sample in the sample vial. In contrast with hydrodynamic injection the sample only needs to touch the capillary tips.

Short electrodes (**G7100-60033**) in combination with hydrodynamic injection are used to minimize sample carry-over to other vials. At appropriate low-fill volumes present in the sample vial during injection, electrodes are not submerged. With this procedure there is no interaction between sample and electrode, such as electrochemical reactions or adhesion of sample to the electrode.

NOTE

This precaution can be programmed into the timetable by a combination of the commands "load vial" and "wait".

Using an Injection Program

The injection table is used for advanced injection tasks such as:

- Injection from different vials, for example ITP.
- Multiple injection modes, for example spiking.
- Injecting a buffer plug after the sample to prevent sample loss after applying voltage.

The online help provides more information on this.

Timetable

In the **CE Timetable**, present on the right-hand side of the CE **Method Setup** dialog you can choose to set and change values for the following parameters at a certain time during the run:

- Voltage
- Current
- Power
- Pressure
- Inlet vial
- Outlet vial
- Cassette temperature
- Lower alarm limit for current
- High pressure (only in CEC mode)

Time programming of these parameters can be useful for example to:

- Define gradients (for voltage or temperature).
- Change the buffer during the analysis (inlet and outlet).
- Apply pressure during the analysis.
- Change the outlet vial during analysis.
- Change the alarm limit for current for example, when changing the buffer system during the analysis.

Instrument Curves

Using a further tab in the CE **Method Setup** dialog box you can choose to monitor and store the following CE-specific raw data. If selected, they are stored in addition to the detector signals defined:

For the CE:

- Leak current
- Internal pressure
- Replenishment level pressure
- Vacuum
- Injection pressure
- External pressure
- Tray temperature
- Cassette temperature

For the DAD:

- Board temperature
- Optical unit temperature
- UV lamp anode voltage

Detector Parameters

The parameters for the detector are set in the DAD tab of the **Method Setup** dialog used in the Agilent ChemStation (see [Figure 20](#) on page 107). On the left hand-side you can define the following settings:

- Used signals, detection and reference wavelengths, bandwidths
- Peak width or detector response time
- Detector stoptime and posttime

Signals, Wavelengths and Bandwidths

Detector parameters allow you to define the detector signals you want to acquire during the analysis and store in the data file. For each signal the wavelengths and bandwidths of sample and reference wavelengths are defined.

It is recommended to do runs without using reference wavelengths to minimize baseline noise. Reference wavelengths can help to compensate for signal drifts. As reference wavelengths are usually measured at higher wavelengths, the baseline noise is increased when using reference wavelengths. Therefore, it might be appropriate to record the two signals at the same wavelength; one using the reference wavelength and one without.

The signal you want to store in the data file is displayed below the lamp icon of the DAD tab of the instrument diagram.

Peak Width

The peak width sets the optimum response time of your detector. If you have similar peak widths in your electropherogram, set it to the width of one of the peaks in your electropherogram. If the peak widths are very different you can use the detector timetable to adjust the peak width setting during the analysis. If your peak width setting is too narrow, peaks might be not detected automatically. This will lead to no spectra for the wider peaks being acquired and in case of peak dependent detection.

Spectra

On the right-hand side of the DAD **Method Setup** dialog an expandable page for advanced settings is visible.

Under **Spectrum** you can define at which points on a signal, spectra will be taken and saved. You can choose to acquire them:

- At the upslope, apex and downslope of the peak
 - All in a peak
 - All spectra during analysis
 - Every 2nd spectrum
- or
- None

You can combine this setting with a wavelength range. By restricting the wavelength range to the range of your interest you can save some disk space; the size of your data files will be smaller because the full spectrum is not stored.

A peak height threshold avoids spectra that are taken from small peaks.

If spectra are recorded, a spectra icon is visible in the detector diagram of the graphical user interface.

Detector Stoptime and Posttime

You can define a special stoptime and posttime for your detector. If you want to use the times set for the Agilent CE instrument, set them to as CE.

Analog Output

The **Analog Output** section enables you to set up the zero offset and attenuation of the analog output signal.

Indirect UV

The section **Indirect UV** enables you to set a margin for peaks with negative absorbance. Set it as close to the baseline as possible but allowing enough range for the negative peaks you expect to get. The higher the value, the lower the resolution of the detector, and the greater the baseline noise. Set this value only if you expect negative absorbances greater than -100 mAU.

Autobalance

Enables you to zero the baseline of the diode-array detector before or after an analysis. You can use the autobalance feature to ensure that the baseline is reset to zero either before the Prerun time has elapsed, or after the Postrun time has elapsed.

Lamps On Required for Acquisition

Deselect the parameter **Lamp on required for analysis** if you also have a different detector configured than the built-in DAD. Such a detector could be an MSD (mass sensitive detector), LIF (laser induced fluorescence) or a CCD (contactless conductivity detector). This option allows you to save lamp life time.

Detector Timetable

Using the **Timetable** for the detector, you can program the time for the following parameters during your analysis:

- Measuring wavelength for signals A to H,
- Measuring bandwidth for signals A to H,
- Reference wavelength for signals A to H,
- Reference bandwidth for signals A to H,
- Spectra acquisition mode,
- Peak width
- Threshold.

By programming the time for these parameters you can optimize the required disk space for your spectra by switching the spectra acquisition on and off. Thus, you can acquire spectra for parts of the electropherogram only.

Switching the Lamp On

The lamp can be switched on automatically as part of initialization (see “DAD Control” on page 200).

To switch the lamp on or off manually:

- 1 In the DAD diagram click right on the lamp icon. A menu will offer the option to switch the lamp on or off depending of the current status. Please see [Figure 23](#) for the different lamps states in the DAD diagram.

Figure 23 UV lamp status Off, Ignition, On

- 2 To ignite the lamp from the menu go to **Instrument > More DAD > UV Lamp** and select **Switch Lamp On** or **Switch Lamp Off**. Please see [Figure 23](#) for the different lamps statuses in the DAD diagram.
- 3 Prior to a run, allow the lamp to warm up for about 1 hour.

If the warm up time of the lamp is too short, a baseline drift in the electropherogram may occur.

Preparing the Replenishment System if Needed

During multiple analyses, the composition of the buffer can change. One way to refresh the buffer is replenishment.

The replenishment system automatically empties and refills the buffer vials in the tray with new buffer from the electrolyte bottle. Typically, replenishment is done before every or after every 3 to 4 runs, depending on the stability and volume of the running buffer.

If you do NOT want to use the replenishment system, you still have to put empty bottles (electrolyte and waste) into the replenishment system at the bottom of the instrument. The bottles are also used as a pressure and vacuum reservoir. Missing bottles lead to pressure and/or vacuum error messages.

WARNING

System under pressure

Pressure is applied to the bottles.

- Only use bottles with part number **9301-6531** or **9301-6532** with this instrument. These bottles are certified to be used under pressure (check the bottles for any scratches or cracks and replace them if necessary). Older instruments were shipped with the bottles **9300-1748** or **5042-6478**. These bottles have a protective coating and can still be used as an alternative.

If you want to use the replenishment system, you have to fill the electrolyte bottle with buffer. Use about 100 mL to make sure the frit is submerged in buffer. Do not fill more than 400 mL buffer into the bottles to ensure the system can work properly. When using buffers containing detergent (for example SDS), the waste bottle should contain some liquid (for example a few mL of isopropanol) to destroy the detergent bubbles. Detailed information on buffers is given in the primer *High Performance Capillary Electrophoresis (5990-3777EN)*, delivered together with the instrument or numerous application notes available on our web page.

Make sure the waste bottle is empty so it can hold the same amount of liquid as in the electrolyte bottle.

When using the smaller 100 ml electrolyte bottle, only use this for the electrolyte. Do not use the 100 ml bottle as the waste bottle.

CAUTION

Instrument damage by corrosion

Components of the replenishment system can be damaged by aggressive or corrosive buffers, organic solvents or evaporated components from these.

- Avoid using aggressive or corrosive buffers or organic solvents that can damage the liquid handling system. Buffers for replenishment or evaporated components from these buffers will come into contact with the following material: polytetrafluoroethylene (PTFE), stainless steel (SST, Grade 316), fluorine rubber (FPM, Viton), polyaryletheretherketone (PEEK), perfluoro-ethylene-propylene (FEP), glass (filter + bottles). A typically acceptable pH range is pH 3-11; for example formic acid can be used in concentrations up to 0.1 %.
 - If you would like to use a buffer composition exceeding the given conditions, we recommend running an analysis without using the replenishment system but individual vials.
-

Before Filling the Replenishment System

The replenishment system (bottles and tubing) should be cleaned when:

- Using it for the first time.
- Changing the replenishment buffer.
- If the replenishment system will be idle for some time.

Filling the Replenishment System

To prepare the replenishment system.

- 1 Click one of the replenishment bottles in the CE diagram.
- 2 Choose **Change Bottles** to release the pressure before you open the bottles.
- 3 Open the front door at the bottom of the front panel.

WARNING

Hazardous organic solvents and sample residues

The waste may contain organic solvents and residue from your sample.

- Wear rubber gloves and safety glasses when handling electrolytes and deposit the waste in a safe waste container.
-

- 4 Take out the electrolyte bottle and unscrew the cap. The electrolyte bottle has a frit and tubing inside. Avoid contaminating the frit (see [Figure 24](#) on page 127).

Filling the Replenishment Bottles

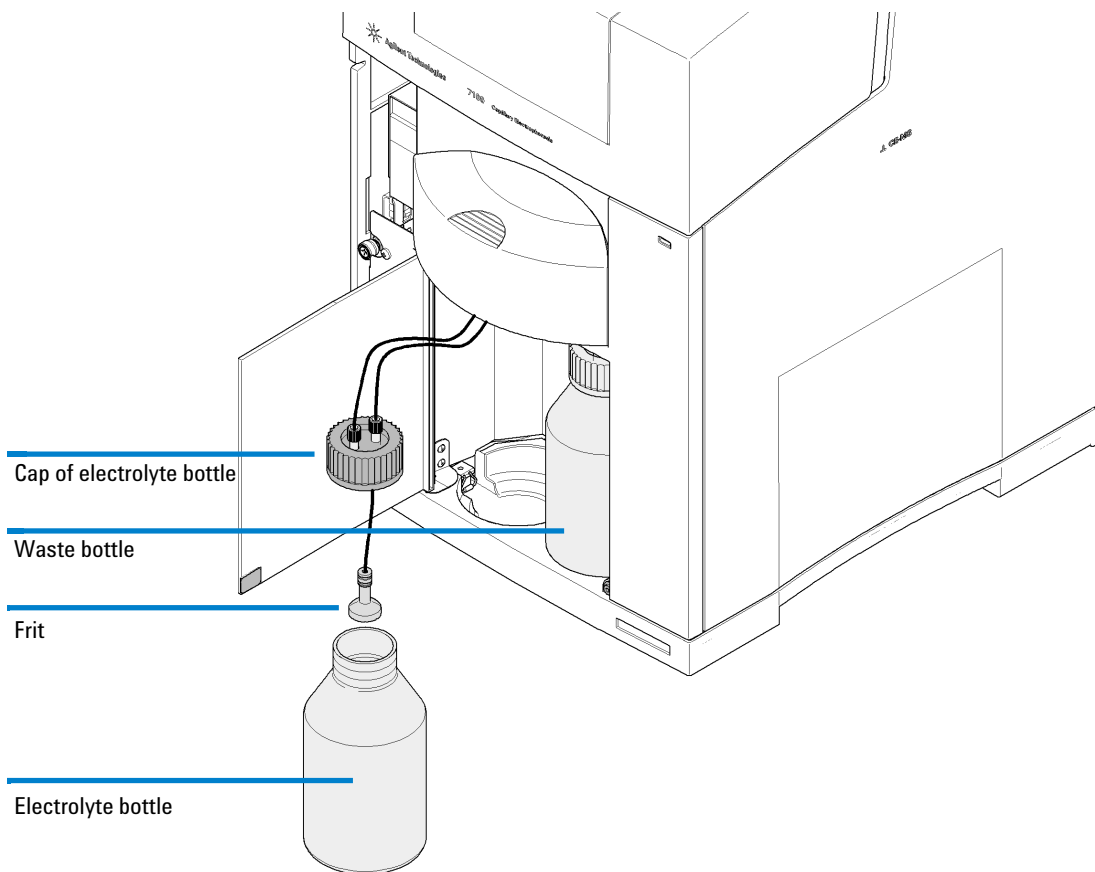
Figure 24 Filling the replenishment bottles

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

1 Make sure the buffer is filtered; use a 0.25 µm filter. For buffers with a high viscosity, a larger pore size, for example, 0.45 µm is also suitable.

2 Fill the electrolyte bottle with buffer.

For 500 mL bottles (part number **9301-6531**), use a minimum of 100 mL to make sure the frit is submerged in buffer. Use a maximum of 400 mL. For 100 mL bottles (part number **9301-6532**), use a minimum of 40 mL to make sure the frit is submerged in buffer. Use a maximum of 80 mL.

3 Tighten the cap.

Make sure the cap is tightened firmly, as pressure will be applied to the bottle.

4 Choose **OK** in the **Change Bottles** dialog box.

Wait until the pressure has built up again, that is the instrument reaches the ready state.

The right-hand bottle is the waste bottle. It is used to store buffer emptied from vials in the tray. Empty the bottle whenever you refill the electrolyte bottle, following the same procedure as described above. Empty the bottle frequently. The error message "Waste Bottle Full" appears if the waste bottle is full.

Preparing Vials for Buffer and Sample

Vial Handling

- Only use only recommended vials (see Table 4) when working with the Agilent CE instrument. Vials must be suitable to be pressurized to avoid busting vials (glass splinter can damage the lifter and block the tray).
- It is recommended cleaning the vials prior use (flush them with water and some alcohol, use compressed air to dry them), otherwise particles may enter the capillary (current problems or blocked capillary).
- All vials must be capped. Use recommended caps only, see [Table 13](#) on page 130.

WARNING





Instrument damage

Avoid damage to the pre-puncher, electrodes and lifter.

- Do not use crimp caps or any other metal caps when working with the Agilent CE instrument.
-
- Make sure there are no air bubbles in the vials, especially when using microvials for samples. Current problems may occur or, in the worst case, the capillary may break (heat generation or no conductance to the outlet buffer).
 - Ensure an adequate liquid level for sample, buffer and waste vials. See [Figure 25](#) on page 131.

5 Using the Agilent 7100 Capillary Electrophoresis System
 Capillary Electrophoresis, CE

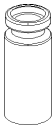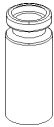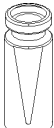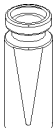Table 13 Recommended vials and caps for the Agilent CE instrument

Item	Description	Part Number	To Be Used For
	Vial clear glass 2 mL Pack of 100	5182-9697	Sample and buffer
	Pack of 500	5183-4623	
	Vial amber glass 2 mL Pack of 100	5183-4619	
	Vial polypropylene 1 mL Pack of 100	5182-0567	Sample and buffer
	Vial polypropylene glass lined 100 µL Pack of 100	9301-0977	Sample only
	Vial polypropylene 100 µL Pack of 1000	9301-0978	Sample only
	Snapcap polyethylene olefin septum for chemical resistance (‘white’ color) Pack of 100	5181-1507 5181-1513	Sample and buffer, but not for multiple use
	Pack of 500		
	PUR polyurethane septum. Preferred standard caps (‘clear’ color) Pack of 100	5181-1512	Sample and buffer

The recommended maximum liquid level in the vials is 1.8 cm measured from the bottom of the vial (see [Figure 25](#)). The minimum level depends on the type of liquid the vial is used for (sample, buffer or waste).

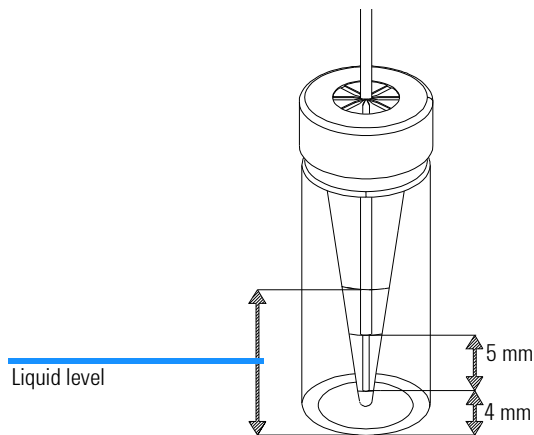
Figure 25 Liquid level and lift offset for standard electrodes

NOTE

To touch the standard electrodes (part number **G7100-60007**), the liquid level must comply with the minimum condition:

Liquid Level = L = Lift offset + 5 mm up to electrode + 1 mm guard band = 1 cm. See [Figure 25](#).

For short electrodes (part number **G7100-60033**), the liquid level must comply with the minimum condition:

Liquid Level = L = Lift offset + 10 mm up to electrode + 1 mm guard band = 1 cm.

To touch the capillary, only the liquid level must comply with the condition for both types of electrodes:

Liquid Level = L = Lift offset + 1 mm guard band = 0.5 cm

Sample Vials

The level of sample in the vials depends on the injection technique used. Please see the section “[Injection](#)” on page 117.

Hydrodynamic Injection

When using injection by pressure, the sample level only needs to touch the capillary. The minimum liquid level in the sample vials for this technique is 5 mm for standard and short electrodes. The short electrode filling height should not exceed 6 mm. This electrode is designed to minimize take-over of sample that absorbs to or simply coats the electrode.

Electrokinetic Injection

When using injection by voltage or current, the sample level needs to touch the electrodes. Therefore, the minimum liquid level in the sample vial is 10 mm. Short electrodes (**G7100-60033**) are not suitable for electrokinetic injection.

NOTE

The effective length in capillary electrophoresis, defined by the distance from the injection end to the detector, generally ranges from 40 to 75 cm. Much shorter effective lengths, such as 8.5 cm from outlet side to the detector are also sometimes very useful. Short effective lengths have the advantage of reduced analysis times, as well as improved peak efficiency and sensitivity.

Hydrodynamic or electrokinetic injection from the short end of the capillary can be achieved by placing the sample vial at the outlet, the buffer vial at the inlet, and then applying a vacuum or a respective electrical field. With the Agilent CE system, up to 100 mbar pressure or vacuum can be applied to the inlet to inject sample at the inlet and outlet, respectively.

Buffer Vials

When the run is started, the liquid level in the vial should touch the electrodes to enable a current to flow. Therefore, the minimum liquid level in the buffer vials is 1 cm for standard electrodes and 1.5 cm for short electrodes. A buffer level of 1.4 cm is recommended for standard electrodes and 1.8 cm for short electrodes. You can use the replenishment system to fill the buffer vials. You can also use the replenishment system to precisely lower the liquid level for all buffer vials to an equal height if you overfill them initially.

Outlet Vials

Outlet vials (for example, waste vials) should also be filled with enough liquid to touch the capillaries. Thus, the liquid flow through the capillary does not produce droplets at the capillary end, which can cause pressure. This would influence the injection amount. If the waste vial is filled with liquid, outgoing droplets cannot be drawn up the electrode by capillary effects but are flushed completely out of the outlet electrode. Therefore, they cannot contaminate the lift head or the next outlet vial. Outlet vials that are used during analysis (for example, for fraction collection) need to be filled to a higher level (1 cm). The liquid must touch the electrode to enable a current to flow.

Vial in Position 49 (Replenishment Lift)

Using replenishment, it is good practice to have a capped vial filled with buffer or water in position 49 (= home position of the replenishment lift). The replenishment needle stays in the liquid and is prevented from buffer precipitation and the resulting blockage of the needle. Water in vial 49 can change the buffer composition inside the replenishment needle and the adjacent tubing over time due to diffusion. This effect might influence the reproducibility of analysis. Therefore, we recommend using buffer.

Inserting the Vials

The tray of the Agilent CE instrument is used for both buffer and samples. The CE diagram of the Agilent ChemStation controls the tray.

- 1 From the CE diagram open the control menu by clicking right at the tray icon.
- 2 The dialogue **Get Vial...** allows you to enter the vial number or tray position.
- 3 Wait until the tray has stopped moving.
- 4 Open the tray door.
- 5 Access the vial which is in the front.

The number is engraved in the sample tray.

NOTE

For safety reasons, the tray can only be moved by the Agilent ChemStation when the tray door is closed.

- 6 Click **Done** after you have accessed all the vial positions you need.

The tray will perform a complete turn every time the tray door is closed in order to read-out the occupied tray positions with an internal vial sensor. Methods will only run if all specified vials are present. The vial sensing procedure might also be delayed if any operation prevents the vial handler temporarily from turning.

Manual Operation

If you want to operate the tray manually or access positions of the tray, it is recommended that you remove the vials from the lifts first. This avoids the need to fill the positions for those vials in the tray again. Use the **Unload** option in the individual lifter context menus to remove the vials from the lifts.

During the run, you can access the vial tray any time the software is not trying to access it. If you try to access the tray during the run, you cannot remove the vials from the inlet and outlet lifts. Therefore, you have to be careful not to fill the positions for those vials in the tray again. Check in the CE diagram which vials are currently located in the lifts.

NOTE

If you insert a vial in a tray position that is already occupied, but where the vial is currently loaded to a lifter, a message will request that you remove the duplicate vial. A duplicate vial is brought to the front position.

Positioning Vials for Sample and Buffer

- 1 Note the positions of the vials in the table (see “[System Vialtable](#)” on page 104).
- 2 Fill the sample into a sample vial and put the vial in the tray, for example in position 7. This position is used in the **Sample Information** screen or the **Sequence Table** screen later.
- 3 Fill at least two vials with the buffer you want to use for your analysis.
- 4 Position those vials in the sample tray. For example, use positions 5 and 6. Write down the positions. When creating the Agilent CE method you must enter the positions of the buffer vials as CE **Home Values** for the inlet home and outlet home vial. After the sample has been injected, those two vials are positioned at the two electrodes.
- 5 When using the replenishment system, you can start off with empty buffer vials and use the replenishment system to fill them automatically with buffer.

Removing the Capillary Cassette from the Agilent CE Instrument

Your Agilent CE instrument comes with an empty capillary cassette installed.

To remove the capillary cassette from the Agilent CE instrument, you have to do the following steps:

- 1 Choose the cassette icon in the CE diagram and select **Change Cassette...** from the menu. This removes the buffer vials from the ends of the capillary and places them back to their position in the tray. This will take approximately 5 seconds.
- 2 Open the top cover to access the cassette. Press a spring located in the middle of the front lid to release the latch. [Figure 26](#), “Position of installed cassette,” on page 135 shows the position of an installed cassette.

NOTE

If you do not lower the vials from the loaded position, buffer, which may be transferred via the capillary into the electrode inserts, precipitates and may block the electrodes or damage the capillary.

- 3 Move the cassette guide to the front.
- 4 Carefully pull the cassette out of the cassette guide. See [Figure 27](#) on page 136.

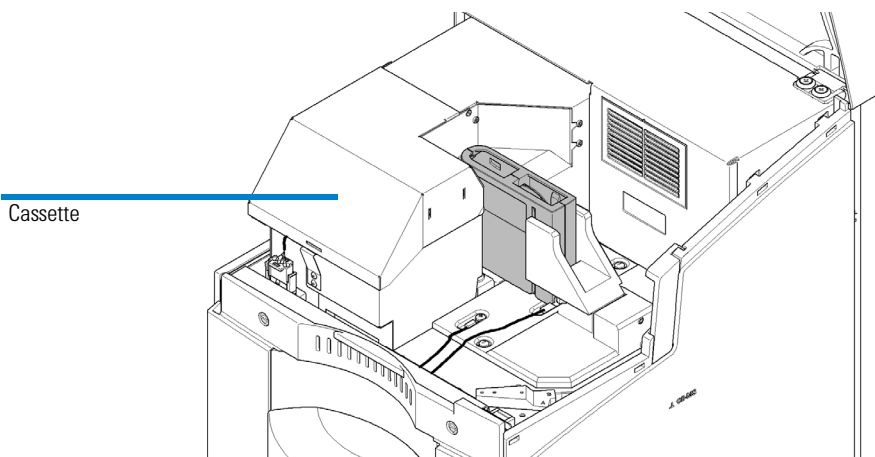
Figure 26 Position of installed cassette

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

NOTE

The DAD detector is equipped with a white LED located near the outlet electrode. This light will assist you when inserting the capillary.

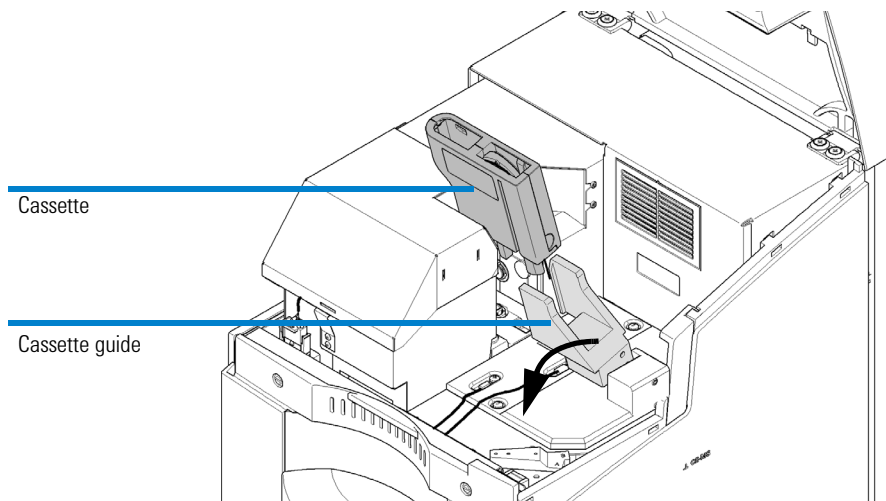
Figure 27 Removing the cassette

Installing a Capillary in the Alignment Interface

The following procedure describes how to install a capillary into an alignment interface. A capillary must have an alignment interface installed before it can be installed into the capillary cassette in order to align the inner capillary with the light path of the diode array detector

Material Needed to Install a Capillary in the Alignment Interface

- Capillary
- Corresponding alignment interface
- Capillary insertion tool (separate item) or part incorporated in cassette
- Safety glasses

WARNING

Risk of injury of sharp capillary ends.

Sharp capillary ends can injure your eyes or skin.

- Wear safety glasses when opening the cassette and handling the capillary to protect your eyes and body.

NOTE

Alignment interfaces for the 7100 CE instrument differ from alignment interfaces for earlier instruments. Interfaces from earlier instruments do not fit into the diode array detector block.

Color Coding of Capillaries and Alignment Interfaces

The alignment interface must correspond to the type of capillary used. There are special alignment interfaces for straight capillaries and for extended light path capillaries. The alignment interface must also correspond to the inner diameter of the capillary. To make it easier to match the capillary and corresponding alignment interface, we use the following color coding on both the sleeve of the alignment interface and the plastic stop on the capillary.

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

Table 14 Alignment interfaces

Sleeve Color	Part Number	Used For*
Black	G7100-60150	25 µm id capillaries with extended light path
Green	G7100-60210	50 µm id straight capillaries
Red	G7100-60230	50 µm id capillaries with extended light path
Blue*	G7100-60310	75 µm id straight capillaries
Yellow	G7100-60330	75 µm id capillaries with extended light path
Green	G7100-60400	50 µm and 75 µm capillaries for CE/MS applications

* Use the blue alignment interface for 100 or 150 µm id capillaries

- 1 Make sure the color of the sleeve in the alignment interface matches the color of the plastic stop on the capillary.
- 2 Remove the protective plastic cap that covers the detection window.

NOTE

The capillary is very fragile at that point and it is important to keep the detection window clean.

- Do not touch the capillary at the detection window or the capillary tips, where the polyimide is burnt off.
- Clean the detection window and the capillary tips as described. Put the insertion tool on a flat surface, for example, your lab bench.

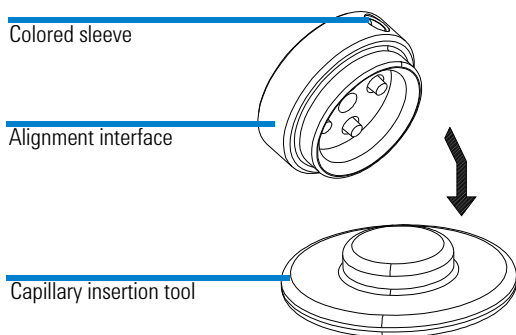
Figure 28 Alignment interface and insertion tool

- 3 Put the alignment interface on the insertion tool and press down firmly. Press only on the outer ring to release the spring that secures the capillary in the alignment interface.

NOTE

Four plastic pins in the alignment interface belong to a feather-driven mechanism to fix the capillary tight in an optimal position for the light path. Press these pins with the alignment tool when inserting a capillary. Insertion attempts of capillaries without this can damage capillary ends and may lead to dissatisfying results.

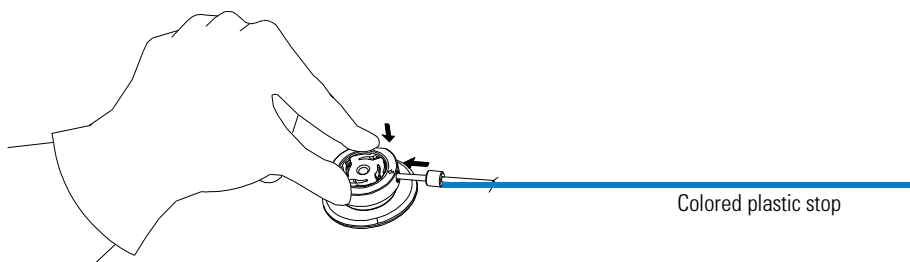
Figure 29 Alignment interface and insertion tool

- 4 Slide the capillary through the opening in the direction of the arrow printed on the interface.
- 5 The capillary has a plastic stop fitted. Pull the capillary through the alignment interface until the stop is reached.

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

- 6 While still slightly pulling the capillary, stop against the alignment interface, and release the alignment interface from the installation tool to secure the capillary in the alignment interface.

The detection window or the extended light path of the capillary must be visible through the detection interface. If aligned correctly, you can see the slit of the interface behind the detection window. See [Figure 30](#) on page 140.

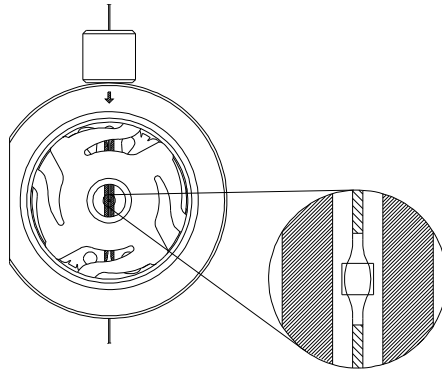
Figure 30 Extended light path aligned correctly

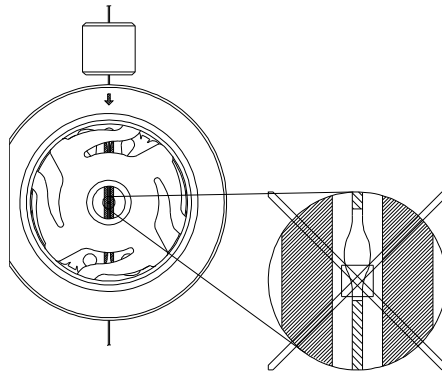
Figure 31 Extended light path aligned incorrectly

Installing a Capillary in an Empty Capillary Cassette

This procedure describes how to install a capillary in an empty capillary cassette. If the cassette has already been installed in the Agilent CE instrument, remove it as described above.

What You Will Do

- 1 Open the capillary cassette.
- 2 Install the capillary in the empty cassette.
- 3 Close the capillary cassette.

WARNING

Risk of injury of sharp capillary ends.

Sharp capillary ends can injure your eyes or skin.

- Wear safety glasses when opening the cassette and handling the capillary to protect your eyes and body.

Open the Capillary Cassette

- 1 Place the cassette flat on a solid surface with stabilizing pins to you.
- 2 Hold the bottom part (for example with both of your thumbs) and slide the top part 8 mm (0.3 inch) gently towards you.
- 3 Lift and remove the top part.

NOTE

Do not use excessive force to open the cassette. This might break parts inside the cassette.

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

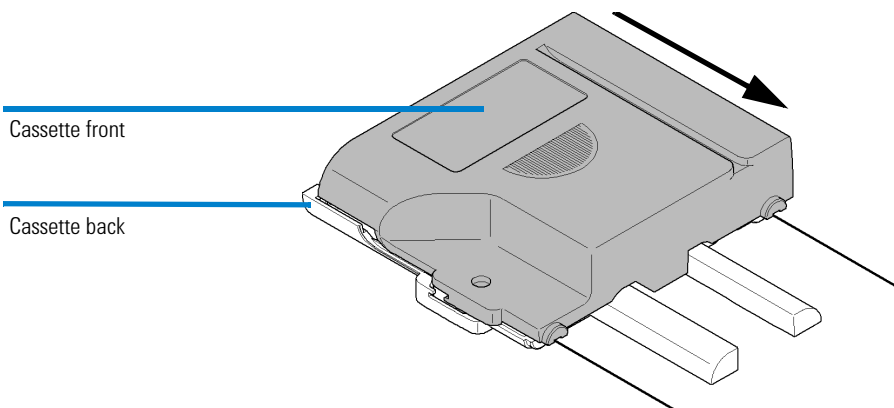
Figure 32 Opening the cassette lock

Installing the Capillary

- 1 Hold the capillary at the alignment interface.
- 2 Insert the alignment interface into the alignment interface holder on the cassette. Make sure the flat side of the alignment interface is at the outer edge of the cassette, see [Figure 33](#) on page 143.
- 3 Make sure the capillary slides into the capillary guiding slit and a guiding hook above the area of the alignment interface.
- 4 Wind the capillary around the reel in the center of the capillary cassette. The pins surrounding the reel have grooves to secure the capillary.

- 5 Start with the inside of the upper pin to prevent the capillary from being squeezed when the cassette is closed.

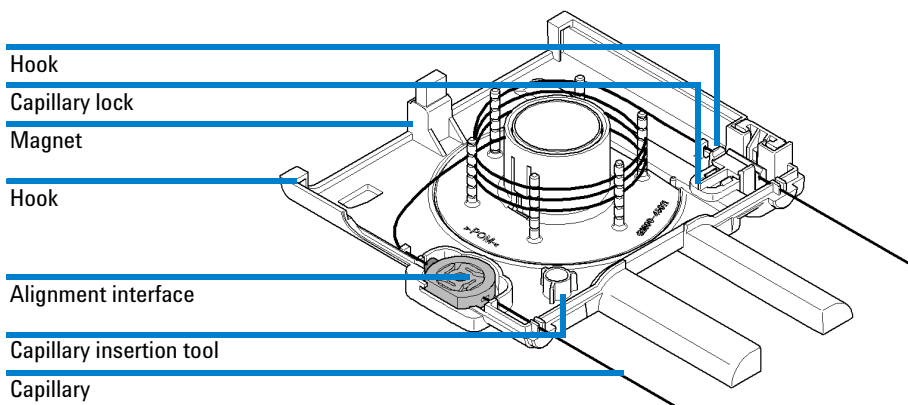
Figure 33 Installing the capillary

Alternate the sides of the pins when inserting the capillary.

NOTE

Make sure the windings of the capillary do not come into contact with each other. As the heat cannot be conducted away at contact points they may get hot and cause separation and detection problems. Such a contact point might also be a predetermined breaking point for arcing.

- 6 Insert the other end of the capillary into the guiding hook on the right.
- 7 Press the white capillary lock towards the center of the cassette and insert the capillary into the capillary lock. Release the capillary lock to secure the capillary.
- 8 Make sure the two capillary ends outside the cassette are the same length as the cassette guiding pins (see [Figure 33](#) on page 143, alignment position) and are parallel to the pins. Use a sheet of paper or another soft item with a straight line to check this. Use the capillary lock to make adjustments on the inlet side.

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

Closing the Capillary Cassette

- 1** Place the top of the cassette onto the bottom part with an offset of 8 mm (0.3 inch).
- 2** Hold the bottom part and slide the top part towards the top of the cassette until it clicks into place.
- 3** Make sure the cassette cover and body form one unit
- 4** Make sure the two capillary ends outside the cassette are both the same length as the cassette guiding pins and are parallel to them. Use the capillary lock to make adjustments to the inlet side of the capillary. The capillary end on the outlet side is fixed; it cannot be moved.

Inserting the Cassette

NOTE

Use the **Change Cassette** task from the CE diagram to access an already installed cassette.

To insert the capillary cassette into the system you have to perform the following steps.

- 1 Choose the cassette icon in the CE diagram and select **Change Cassette...** from the menu.
This removes the buffer vials from the ends of the capillaries and back to their position in the tray. This will take approximately 5 seconds.
- 2 Open the top cover to access the cassette.
- 3 Press a spring located in the middle of the front lid to release the latch.
[Figure 34](#) shows the position of an installed cassette.

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

- 4 Move the cassette guide to the front.

NOTE

Extracting the cassette from the instrument without lowering the lifts will lead to a higher contamination of the inner parts of the electrode or pre-puncher. Crystallization or liquid residue of this contaminations might bias future runs.

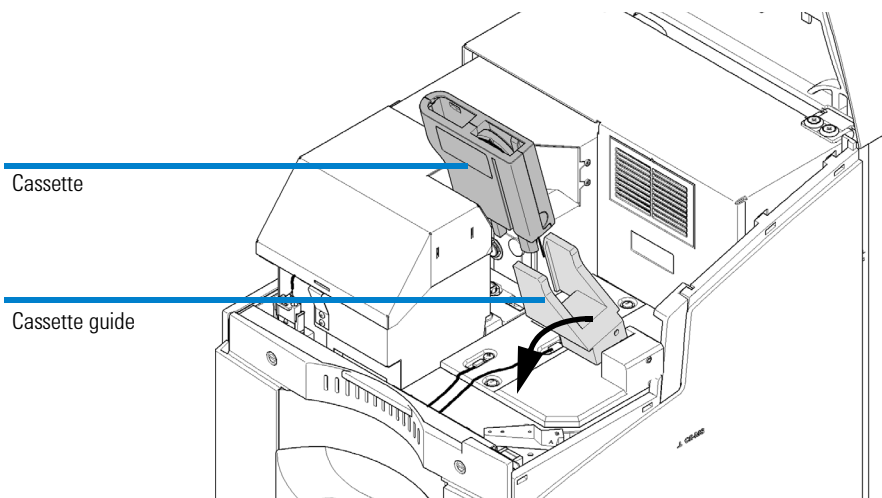
Figure 34 Insertion of cassette

CAUTION

Capillary damage by incorrect insertion into capillary seats

The capillaries should slide directly into the capillary seats. Otherwise, they might break or be damaged.

- Make sure the ends of the capillary are not bent when lowering the capillary cassette.
- If you have problems inserting the capillary into the electrodes, check if they are blocked with salt crystals (see [“Cleaning the Electrodes, Pre-punchers and Insulation Plate”](#) on page 216).

- 5 Slowly lower the cassette until it reaches the insulation plate. While lowering the cassette, insert the capillary ends into the capillary seats of the electrodes.

- 6 Move the cassette guide back again while holding the cassette down. The alignment interface aligns the capillary automatically with the detector.

NOTE

Only alignment interfaces dedicated to the 7100 CE instrument will fit. Interfaces from different instruments will not fit and will lead to unsatisfactory results.

CAUTION

Damage to the cassette

Closing the top cover with an improperly installed cassette may damage the cassette.

- Do not close the top cover, if the cassette is not properly installed.

- 7 Close the top cover and secure it by latching it onto the body of the case. For safety reasons, no power is applied if the top cover is open. Because the cassette was changed, the Agilent ChemStation automatically performs the wavelength calibration in the background. It tests the wavelength calibration of your detector and advises you to make adjustments, if necessary.

If the lamp is off or the cassette is not inserted properly, the wavelength cannot be adjusted correctly.

- 8 Close the **Change Cassette** dialog box by choosing **OK**, enter the capillary information and proceed to setting up a method.

One of the following not ready conditions may appear in the status display of the Agilent ChemStation. Move the mouse over the status display to read the not ready conditions:

- Recalibration necessary
- UV lamp not ready

To correct for these not ready conditions please reinsert the cassette. Make sure the alignment interface snaps onto the ceramic interface of the detector and is not occluded. Make sure the lamp is ignited.

Fraction Collection

In the bioscience and pharmaceutical sectors, the capability of fraction collection is frequently required in order to characterize separated components by off-line analysis techniques such as MALDI/TOF-MS or protein sequencing. CE is in essence a miniaturized separation technique. However there is a continued interest in its preparative possibilities due to its unsurpassed separation efficiency and the fact that it can be performed in free solution.

Bulk production of desired compounds is marginal but the technique remains still attractive for the production of small amounts of purified components which are not available by other separation techniques.

The separated, tiny fractions might be collected for the following purposes:
Sequencing of DNA fragments or peptides,

- Off-line identification by MS and molecular weight determination by MALDI-TOF,
- Establishing biological activity in a single substance, or
- Checking the purity of a collected peak by reinjection
- Small scale preparation of reference substances

The schematic of a sample collection, during a preparative CE run, is illustrated in [Figure 35](#) (5990-3777EN CE primer).

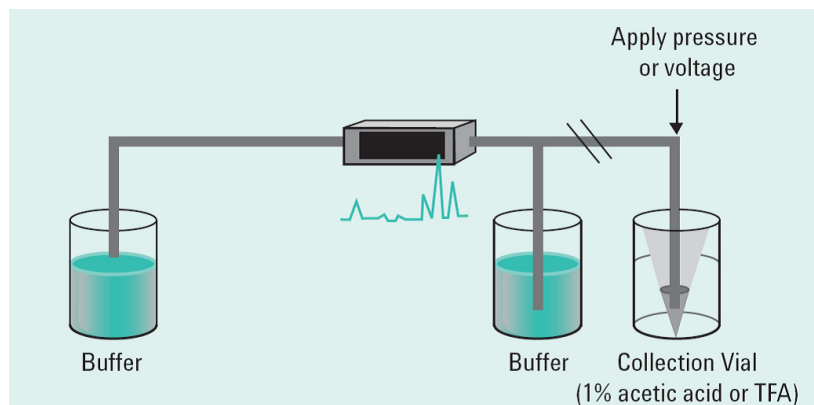
Figure 35 Schematic of sample collections

Fraction collection execution on the 7100 Capillary electrophoresis instrument uses information provided by a built in peak detector. This peak detector resides on the instrument as part of the firmware. From data recorded by the built-in diode array detector the exact time when a peak appears at the outlet of the capillary is calculated. This peak is collected to a collection vial with a window of certain range around it.

Peak detection triggers fraction collection if the signal exceeds certain threshold or a certain slope or a combination of both. Either all peaks in a time window are collected or, for up to four expected peaks, anticipated migration times can be set. The signal of wavelength A is used exclusively.

The length of the capillary between the detection window and the outlet is fixed to 8.5 cm. An algorithm, which is part of the data analysis software (ChemStation C.01.07 SR1 or higher), calculates based on the peak migration time, the elution mode, the capillary length and the electrical field the time in which the peak will approach the end of the capillary. Shortly before this point in time the instrument stops electrophoresis and swaps the outlet home vial to a collection vial. Subsequently either the peak is moved out pressure assisted or by means of electrophoresis for a time defined by the collection width. After this collection step the instrument swaps back to the initial outlet home vial and continues electrophoresis until next fraction collection steps happen or the run stop time elapses.

Major prerequisites for successful fraction collection on the 7100 Capillary electrophoresis instrument is a development of a suitable method that provides stable migration times and peak width and heights of the respective components. Potential applications include fraction collection from capillary zone electrophoresis CZE, micellar electrokinetic capillary chromatography MECC, from capillary gel electrophoresis CGE and even from capillary isoelectric focusing CIEF separations.

Peaks may be collected either by using pressure or electrokinetic elution. Please bear in mind injected amounts are small and peak volumes are generally in the size of nanoliter (nL) volumes. Hence peaks are typically collected into microvials containing minimal volumes of an appropriate solution to have a minimal dilution. Identical peaks are often collected numerous times into the same collection vial. Fill level of the collection vial depends on the collection mode. About 20 μ l liquid are sufficient for pressure elution as only the capillary end needs to be submerged. For electrokinetic elution 50 μ l are required as the outlet electrode needs to be submerged for electrical contact.

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

NOTE

When using short electrodes (part number **G7100-60033**) the outlet vial which is used during fraction collection needs to be filled to a higher level (1 cm). The liquid must touch the electrode to enable a current to flow for electrokinetic elution. Fill level for pressure elution is not affected.

A method suitable for fraction collection employs often a capillary with bigger inner diameter such as 75 μm , 100 μm or 150 μm to maximize sample introduction at injection. This partially sacrifices often resolution or lowers the achieved plate number of a similar analytical run. Goal is to limit heat creation and or time consumption for long runs balanced while achieving separation that is just sufficient for fraction collection. As a consequence buffer composition, voltage, injection, cassette temperature and capillary length need to be optimized differently for preparative tasks.

NOTE

The fraction collection algorithm modifies the CE time table during the run based on actual migration times of peaks during the run. When using fraction collection all timetable entries which use voltage, outlet vial or pressure settings will be deleted from the timetable.

Provide a sufficient number of vials filled with appropriate liquid for pressure or electrokinetic collection mode in consecutive location in the sample tray. The number of vials depends on the number of peaks that are expected to be collected.

The instrument logbook shows the steps done during fraction collection. It provides information when a peak has been detected and if it falls into a search window. The vial into which the peak has been collected and time frame of the collection is noted.

Prior to any fraction collection run capillary information must be provided in the analysis software under **Instrument > Capillaries** (Figure 36 on page 151). In this dialogue capillary information such as the material description and the inner diameter of the capillary is given. The values for total capillary length and effective capillary length, which is the distance from inlet to the detection window, set here are taken into account by the fraction collection algorithm. They are used to calculate migration speed and speed of movements by pressure application and are therefore crucial for successful fraction collection.

#	Installed	Description	Col. Serial#	Batch#	Product#	# Injections	Max. p [bar]	Max. T [°C]	Max. pH	Min. pH	Length	Diameter	Cell	eff. Length
1	YES	Barefused Silica	autoID-40		G1600-60211	30	0	0.0	2.0	0.0	48.5	50.0	Capillary	40.0
2	no	Barefused Silica	autoID-39		G1600-60132	0	0	0.0	2.0	0.0	48.5	25.0	Bubble	40.0
3	no	Barefused Silica	autoID-41		G1600-60232	0	0	0.0	2.0	0.0	48.5	50.0	Bubble	40.0
4	no	Barefused Silica	autoID-42		G1600-60233	0	0	0.0	2.0	0.0	43.5	50.0	Bubble	35.0

Figure 36 Capillary information screen

The Fraction Collection Screen

The Fraction Collection Screen is available from the instruments user interface of ChemStation C.01.07 SR1 or higher. (Figure 37):

Instrument > More CE > Set up Fraction Collection...

It will also automatically appear in the sequence of dialogs for editing entire methods:

Method > Edit entire method ...

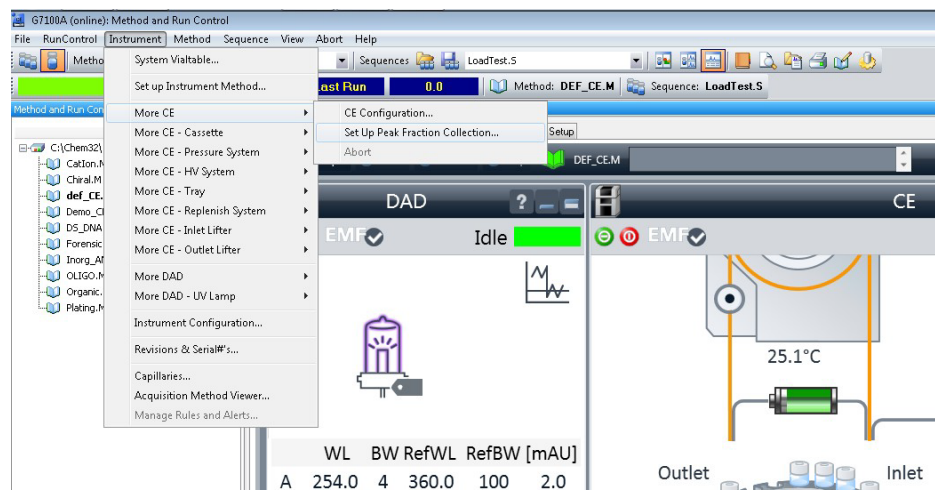
Figure 37 Set Up Peak Fraction collection

5 Using the Agilent 7100 Capillary Electrophoresis System Capillary Electrophoresis, CE

The Fraction Collection Screen is given in [Figure 38](#). The default setting for the Collection mode is OFF. For methods with this setting no fraction collection is done, any value entered in the fields below will be ignored. The content of the dialog depends on the selected Collection mode.

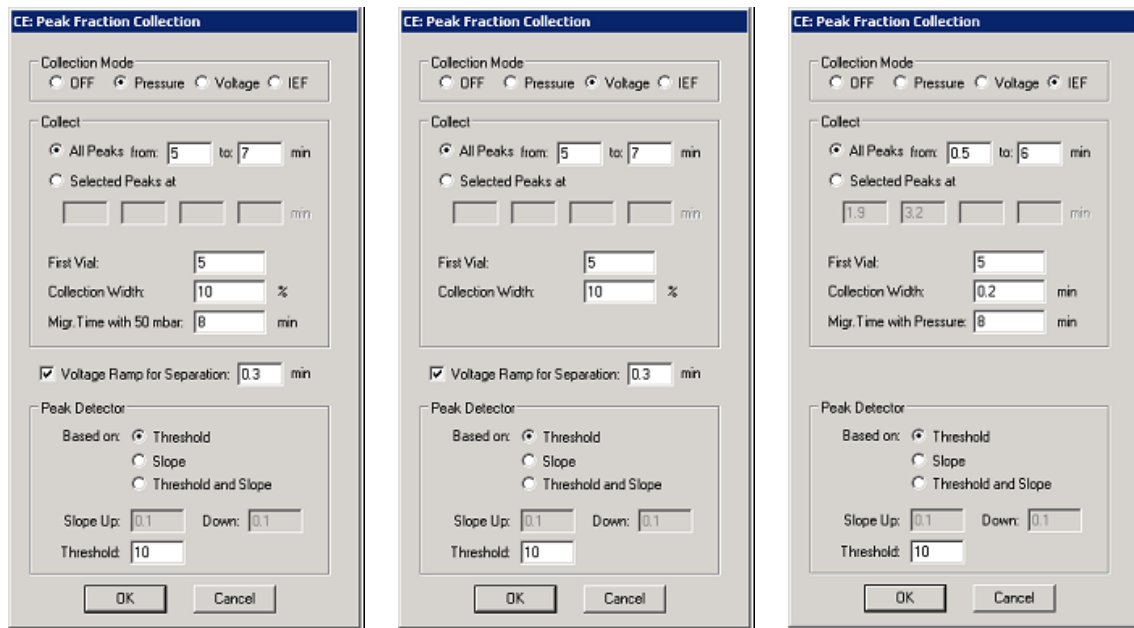
Figure 38 Fraction Collection Screen for pressure (left), elektrokinetic (middle), combined (right) elution of peaks

Collection Modes

Four alternative collection modes are available.

OFF	No fraction collection set up.
Pressure	Moves the peak into the collection vial by applying 50 mbar pressure. The collection vial should be filled with sufficient run buffer or other electrolyte to ensure the leaving liquid from the capillary end is collated with it.
Voltage	Moves the peak into the collection vial by applying the voltage. The collection vial should be filled with sufficient run buffer or other electrolyte to ensure conductance.
IEF	50 mbar pressure is used to move the peak into the collection vial. In the IEF mode the peaks have no electrophoretic mobility so the algorithm to calculate the time when the peaks reach the outlet is different than in collection with pressure only.

Collect

Two alternative options for activation of the peak collection algorithm are available. Depending on the collection mode selected the items in the Collect panel Change.

Option 1	All peaks: from start time to end time in minute
Option 2	Selected Peaks at: mt of four peaks can be set

Select Option 1 (**All Peaks**) if you want to have all detected peaks above a certain threshold collected.

Select Option 2 (**Selected Peaks**) and provide migration times into the fields to collect up to 4 selected peaks expected within each time window specified by the migration time and the Collection width.

For both options enter in the field **First Vial** the start vial where the first fraction is collected can be entered. The second fraction will be collected in the next vial. An error message will be logged in the logbook if the actual vial number exceeds vial position 48. No further peaks will be collected.

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

NOTE

For fraction collection on Selected Peaks the vial handling and elution process consumes time and due to this break in the separation the second, 3rd and 4th peak will migrate at a later time than in runs without fraction collection. In order to collect multiple peaks method development is an iterative process.

NOTE

Do not use the Lower Alarm Limit or time program it when changing the inlet or outlet vial during the run during fraction collection. If Lower Alarm Limit is active the system will stop when changing the vial because the lower alarm limit is reached as the current is zero at this point in time.

Collection Width

The collection width defines the peak collection time window around the peak apex. Depending on the collection mode, it is specified in percentage of the time the peak appears or an absolute time. This is the migration time plus the time the peak needs from the detector to the outlet vial. This is because slower peaks will become broader and therefore need a wider migration time window.

Example: If a collection width of 5% is used for a peak which has an expected migration time of 1 minutes (detection point to capillary outlet-end) then the collected peak width will be 0.6 minutes. Therefore peak collection will begin at 11.7 minutes and continue until 12.3 minutes. If peaks are being collected from an IEF separation, the peak width is not dependent upon migration time so the peak collection time can be defined and remains the same for all collected peaks.

Migration Time with 50 mbar/Pressure

If you use collection by Pressure or IEF the instrument needs to know at which velocity the peaks are moving. In other words, how long it would take for an injected sample plug to move from the capillary inlet to the detector if moved by pressure only. To find this value, create the following method:

- 1 Switch High Voltage off.
- 2 Keep cassette temperature as in the later experiments.
- 3 In the timetable enter: At time 0 minutes internal pressure 50 mbar.
- 4 Set a suitable stop time, for first runs chose of “no limit”.
- 5 Inject your sample or any absorbing liquid (e.g. Caffeine, NaOH, isopropyl alcohol).

- 6 Set a wavelength where your sample or the liquid absorbs (e.g. 235 nm for Caffeine).
- 7 In method step **Preconditioning** flush the capillary with your run buffer for 5 minutes.
- 8 Start the method, read out at what time the sample appears in the detector window.

The time should be entered as "Migr. Time with 50 mbar" or under "Migr. Time With Pressure" for IEF collection mode.

Voltage Ramp for Separation

Sometimes a method starts with a voltage ramp to prevent rapid heating of an injected sample or in order to protect the hardware from damage. The default ramp up rate of the 7100 Capillary electrophoresis is 5 kV/s. Conditions other than this are set in the **Set up CE Timetable** screen. Set a voltage ramp if desired in the section **Voltage Ramp for Separation** of the **Peak Fraction Collection** dialog.

NOTE

The fraction collection algorithm modifies the CE time table during the run based on actual migration times of peaks during the run. When using fraction collection all timetable entries which use voltage, outlet vial or pressure settings will be deleted from the timetable. Therefore set the voltage you want to use in the electric screen and the time for the voltage ramp in the fraction collection screen. The voltage ramp must be completed before the first peak to be collected is expected.

Summary of Steps for Fraction Collection

- Optimize separation method for high peak concentration and migration time reproducibility.
- Fill micro vials with low volume of appropriate solution (see [Table 15](#) on page 156).
- Place micro vials in sample tray (random choice, consecutive order).
- Enter the correct capillary information in the 'capillary information screen.'
- Find and enter the migr. time with pressure values.
- Set threshold or slope according to your peak.
- Run fraction collection method.
- Track detection and fractionation process in online plot and logbook.

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, CE

- Improve method for multiple peaks stepwise.
- Look into the logbook to confirm the fractionation collection steps happened as expected.
- Mix the collected fraction thoroughly when you want to re-analyze the collected fractions.

Be aware that the result depends on many factors like sample concentration, injection time and pressure, capillary length and internal diameter. The online plot of ChemStation allows convenient monitoring the pressure, current and voltage readings of a run. It is highly recommended to track this as well as the log book of the run during method development.

Recommended Collection Modes

Table 15 gives an overview of the recommended collection modes using the various CE separation techniques and proper fraction collection solutions.

Table 15 Collection Modes

Separation Technique	Elution Mode Pressure	Elution Mode Voltage	Collection Liquid
CZE	Preferred	Yes	2% acetic acid
ITP	Preferred	Yes	2% acetic acid
CGE (using gels)	Don't use!	Yes	Electrolyte
MEKC	Yes	Yes	Electrolyte
IEF	Yes	No	Carrier ampholyte

Capillary Electrophoresis, Plus High Pressure CE+p

The Agilent CE instrument is capable of accepting an external gas pressure in addition to the internal pressure. The external high pressure source can be up to 15 bar (218 psi), and is regulated within the instrument via method settings or with direct instructions from the CE diagram between 2-12 bar. This pressure can be applied to one or both vials at the capillary ends. The high pressure option may be required for capillary isoelectric focusing (cIEF) or if using highly viscous replaceable gels in the CGE.

Requirements

- External pressure supply (< 15 bar) of oil-free air or nitrogen.
- Push-fit connector and gas line (supplied in startup kit).
- Outlet lift connection to the tubing system

NOTE

For external pressure supply, a male adapter fitting the instrument's female adapter and PTFE tubing (1/8" OD x 1/16" ID PTFE tubing) are part of the accessory kit of the instrument. Please add a suitable connector to the external pressure supply.

The inlet for the external high pressure gas is situated at the back of the instrument:

5 Using the Agilent 7100 Capillary Electrophoresis System
Capillary Electrophoresis, Plus High Pressure CE+p

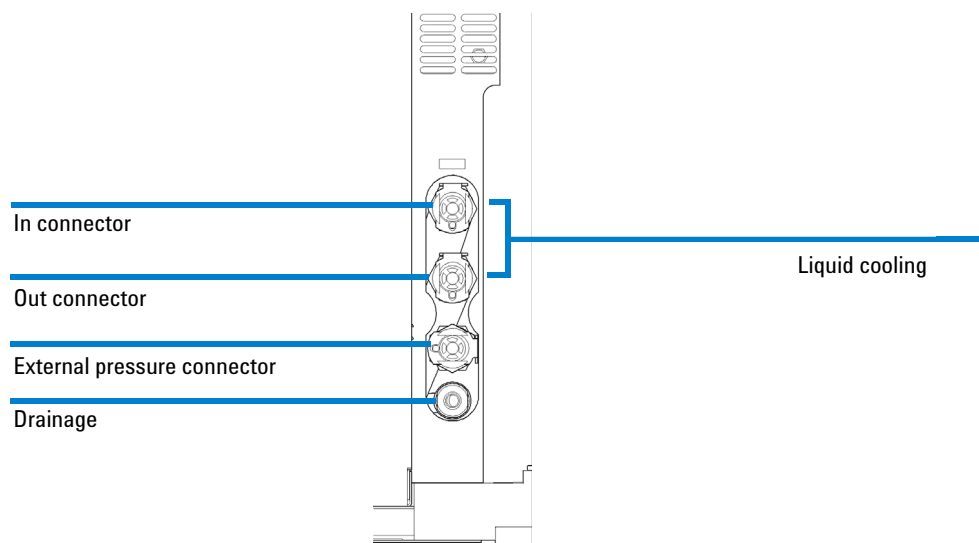
Figure 39 External high pressure gas inlet

Outlet Lift Connection to the Tubing System

The outlet lift is not connected to the tubing and valve system of the instrument by default. Please connect the respective PTFE tube provided in your accessories kit. The tube is labeled with an "8" and has one open end and one fitting.

To check whether this tubing is already installed, you should load a vial in the outlet lifter from the CE diagram and locate the tube coming from the outlet lift. [Figure 40](#) shows an outlet lift with mounted tubing.

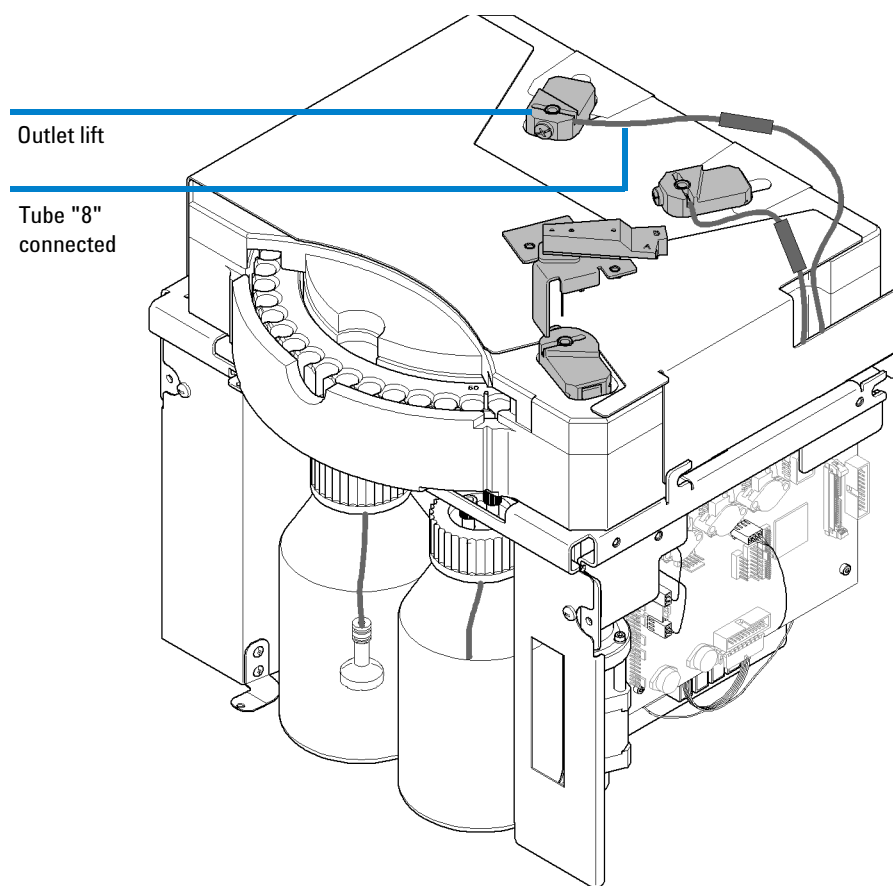
Figure 40 Outlet lift with mounted tubing for high pressure application

WARNING

Damage or contamination of pressure valves.

Make sure tube "8" is connected properly before you apply high pressure to any lift.

High pressure may flush liquid into the pressure valves and lead to permanent damage.

- Connect tube "8" prior to using high pressure.
-

Steps to Install the Outlet Lift Connection to the Tubing

Please refer to the section [“Cleaning the Electrodes, Pre-punchers and Insulation Plate”](#) on page 216 for more details on individual steps of the following procedure.

- 1 Unload all lifter positions.
- 2 Switch off the instrument.
- 3 Disconnect the power.
- 4 Open the front cover by releasing two screws and pulling it towards you.
- 5 Loose two screws to release the liquid handling module.
- 6 Pull the liquid handling module towards you.
- 7 Remove the plastic screw of the outlet lifter (left).
- 8 Extract the pre-puncher from the lift head by moving it straight up with tweezers.

WARNING

Injury by sharp metal edges

Pre-punchers are sharp and might result in injury.

To prevent personal injury, be careful when getting in contact with sharp metal areas.

- 9 Slide the open end of the PTFE tubing over the outlet pre-puncher using forceps or by grabbing it with sandpaper wrapped around it.
- 10 Insert the outlet pre-puncher in the lift head in a vertical movement.
- 11 Tighten the plastic screw.
- 12 Slide the liquid handling module back into the instrument until it is positioned by internal pins.
- 13 Tighten the two screws to fix the liquid handling module.
- 14 Mount the front cover and tighten the two screws to fix.

Preparing the Agilent CE Instrument for High Pressure Use

- 1 Ensure that the gas source is switched off prior to connection.
- 2 Connect the push-fit connector and gas line to the receptor on the rear of the Agilent CE instrument.
- 3 Connect the gas line to the source.
- 4 Switch on the gas supply.
- 5 From the instrument control screen select **Instrument > More CE > CE Configuration**, see [Figure 41](#).

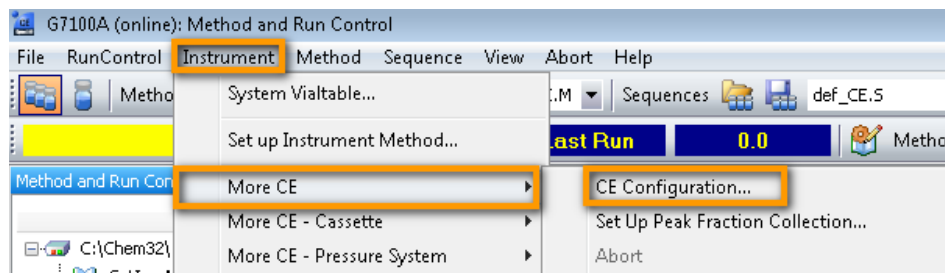
Figure 41 Configuration menu

5 Using the Agilent 7100 Capillary Electrophoresis System
Capillary Electrophoresis, Plus High Pressure CE+p

- 6** Enable CE+p mode by selecting the option **External Pressure Installed**.

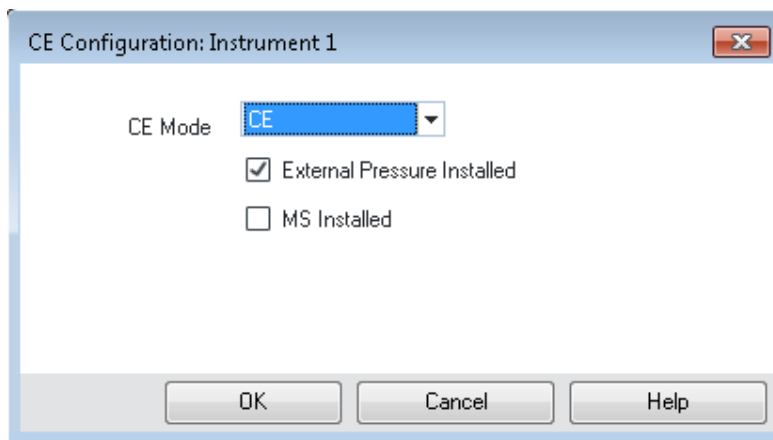
Figure 42 Select CE mode

- 7** Confirm your selection by clicking **OK**.
- 8** Make sure tube “8” is connected, see [“Outlet Lift Connection to the Tubing System”](#) on page 159.
- 9** Once the external pressure is configured a new icon for an external pressure reservoir will appear, see [Figure 43](#).

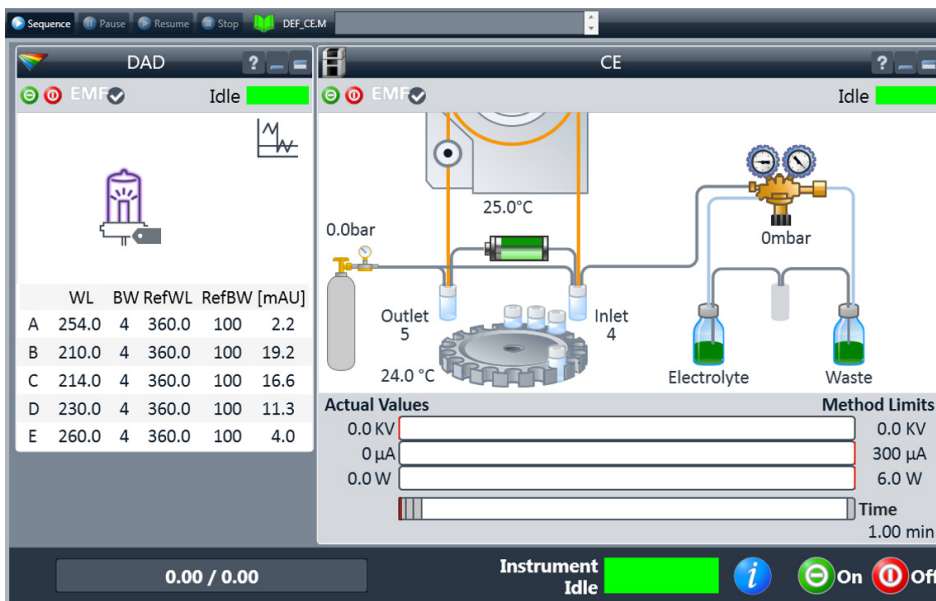
Figure 43 CE diagram with CE+p mode activated

Operation and Function in CE+p Mode

By selecting CE+p mode, some additional functions are available for control via the CE diagram, which will assist in the operation of capillary gel electrophoresis (CGE) with highly viscous buffers. This includes the ability to apply high pressure during preconditioning and postconditioning and the ability to use high pressure to introduce gel or samples to the capillary.

Preconditioning and Postconditioning

When the instrument is operating in CE+p mode, the user can chose to use the **High Pressure Flush** command menu during preconditioning or postconditioning (Figure 45). This means that high pressure will be applied to the capillary inlet vial which may be:

- A set value (2-12 bar)
- Raw external pressure (up to 15 bar): Allows pressurizing the inlet vial and/or the outlet vial. The external pressure up to 15 bar can be applied directly without reduction.

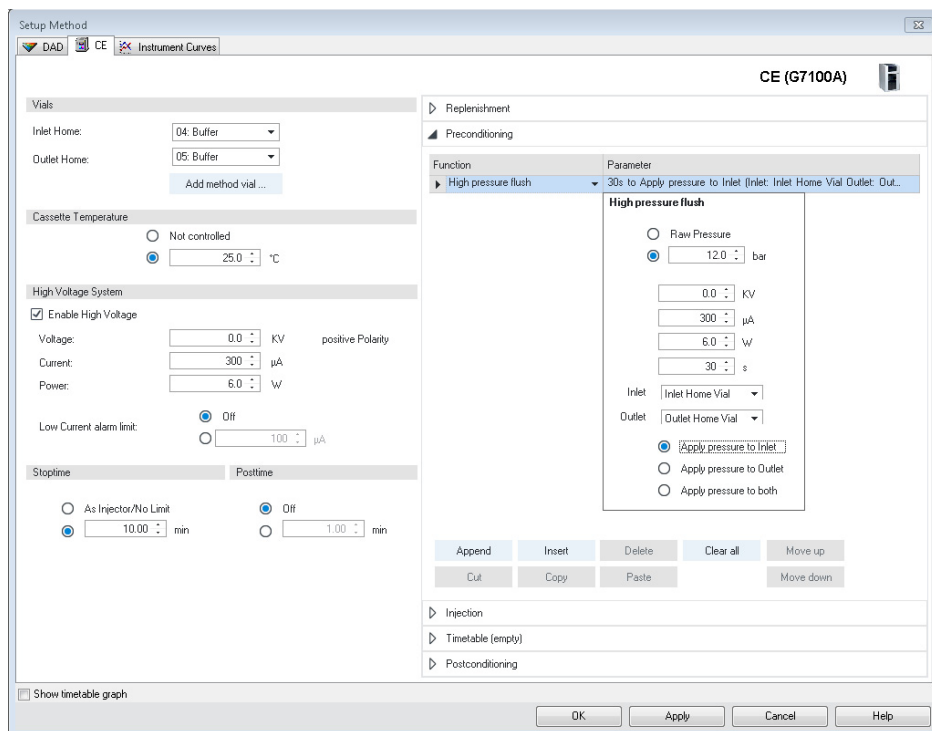
Figure 44 Method setup with CE+p mode activated

5 Using the Agilent 7100 Capillary Electrophoresis System Capillary Electrophoresis, Plus High Pressure CE+p

High pressure flush

Raw Pressure

bar

kV

μ A

W

s

Inlet

Outlet

Apply pressure to Inlet

Apply pressure to Outlet

Apply pressure to both

Figure 45 Parameters for high pressure flush

Other parameters which can be set in the **High pressure flush** command menu include:

- Applied voltage (kV), current (μ A) or power (W)
- Duration (s),
- Inlet vial and/or outlet vial

Injection Options with CE+p

The installed high pressure supply may also be used for injection purposes using the **Timetable** option in the **Injection** field of a method setup.

From this menu you may select the **High pressure flush** command menu during injection with the same parameter choices that are available during Preconditioning.

This may be used to perform a pressure injection while using a capillary filled with replaceable gel.

NOTE

Please note that the external pressure is not as precise as the injection pressure of up to 100 mbar!

Timetable Options with CE+p

The installed high pressure supply may also be used in the general timetable of a method setup. The **Pressure** command is extended by the **External** pressure functionality, see [Figure 46](#) on page 167.

Pressure

- Off
- Internal mbar
- Flush (950 mbar internal)
- Raw External Pressure
- External bar
 - Apply pressure to Inlet
 - Apply pressure to Outlet
 - Apply pressure to both

Figure 46 Pressure options in CE+p mode

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrophoresis, Plus High Pressure CE+p

Timetable controlled changes in the pressure settings are:

- OFF: Pressure usage is switched off.
- Internal, pressure (-100 mbar to +100 mbar): Can be applied to the inlet vial.
- Flush: Standard flush procedure is applied with an internally generated pressure of approximately 950 mbar.
- External: Allows pressurizing the inlet vial and/or the outlet vial. External pressure is reduced to the set value between 2 and 12 bar and applied in a controlled manner.
- Raw External Pressure: Allows pressurizing the inlet vial, and/or the outlet vial. The external pressure up to 15 bar can be applied directly without reduction.

CE+p Control Using the CE Diagram

High pressure may be applied independently of the method for capillary conditioning or cleaning by clicking right at the external pressure icon in the CE diagram (see [Figure 43](#) on page 163). This will cause a parameter setting field to appear (see [Figure 45](#) on page 166). From this field you may choose to apply pressure to the inlet, outlet or both vials, and also whether to simultaneously apply a voltage of -30 to 30 kV for a given time.

Capillary Electrochromatography, CEC

Capillary electrochromatography (CEC) is a fusion of capillary electrophoresis and capillary chromatography. In CEC, solvent is mobilized through a packed capillary column by electroosmotic flow instead of the hydraulic flow used in liquid chromatography LC. Higher efficiencies are achievable because of the uniform flow velocity profile and the ability to use smaller particle sizes in longer columns than is possible with LC. Operation of CEC requires an equal application of high pressure to the inlet and outlet vials in order to suppress bubble formation while avoiding hydrodynamic flow due to a pressure differential.

How to Use Your Agilent CE Instrument for Capillary Electrochromatography

Requirements

- External pressure supply (< 15 bar) of oil-free air or nitrogen.
- Push-fit connector and gas line (supplied in startup kit).
- Outlet lift connection to the tubing system

NOTE

For external pressure supply, a male adapter fitting the instrument's female adapter and PTFE tubing (1/8" OD x 1/16" ID PTFE tubing) are part in the accessory kit of the instrument. Please add a suitable connector to the external pressure supply.

The inlet for the external high pressure gas is situated at the back of the instrument.

5 Using the Agilent 7100 Capillary Electrophoresis System
Capillary Electrochromatography, CEC

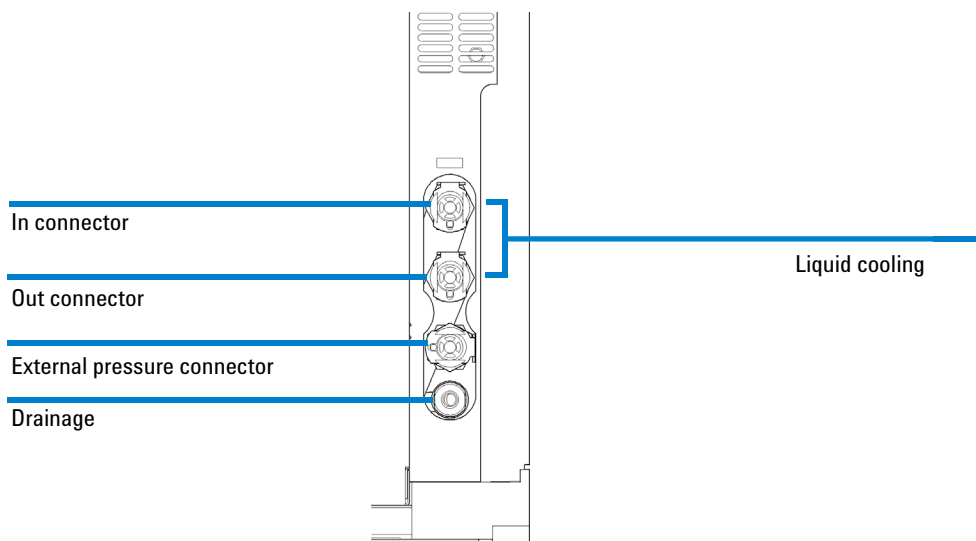
Figure 47 External high pressure gas inlet

Preparing the Agilent CE Instrument for High Pressure Use

- 1 Ensure that the gas source is switched off prior to connection.
- 2 Connect the push-fit connector and gas line to the receptor on the rear of the Agilent CE instrument.
- 3 Connect the gas line to the source.
- 4 Switch on the gas supply.
- 5 From the instrument control screen select **Instrument > More CE > CE Configuration**, see [Figure 48](#).

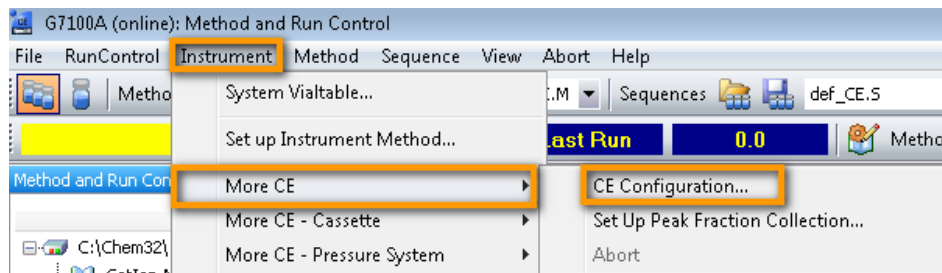
Figure 48 Configuration menu

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrochromatography, CEC

- 6 Enable the CEC mode by selecting the CE mode **CEC** (note: **External Pressure Installed** option is automatically selected and the option is greyed out)

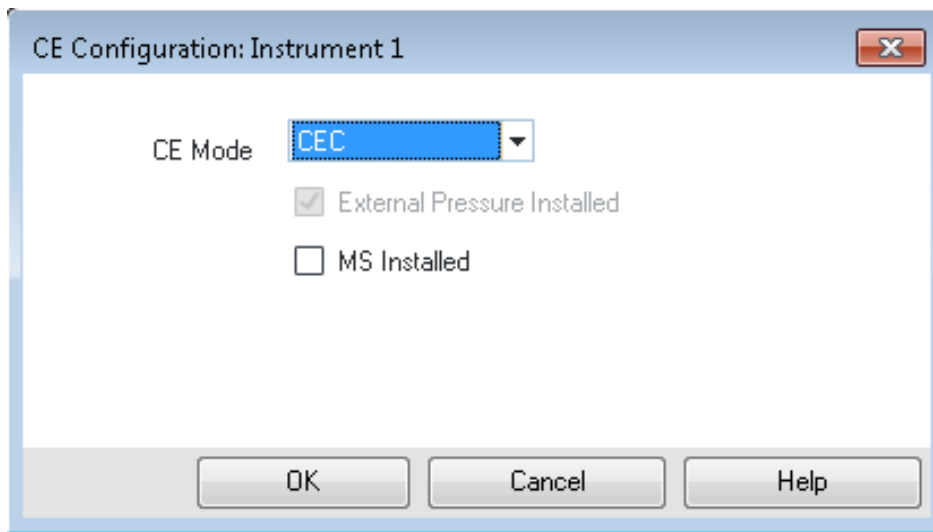
Figure 49 Select CEC mode (to decide which access to config is used)

- 7 Confirm your selection by clicking **OK**.
- 8 Make sure tube "8" is connected, see [“Outlet Lift Connection to the Tubing System”](#) on page 159.

WARNING

Damage or contamination of pressure valves.

Make sure tube "8" is connected properly before you apply high pressure to any lift. High pressure may flush liquid into the pressure valves and lead to permanent damage.

- Connect tube "8" prior to using high pressure.

- 9 Once the CEC mode is configured a new icon for an external pressure reservoir will appear, see [Figure 50](#).

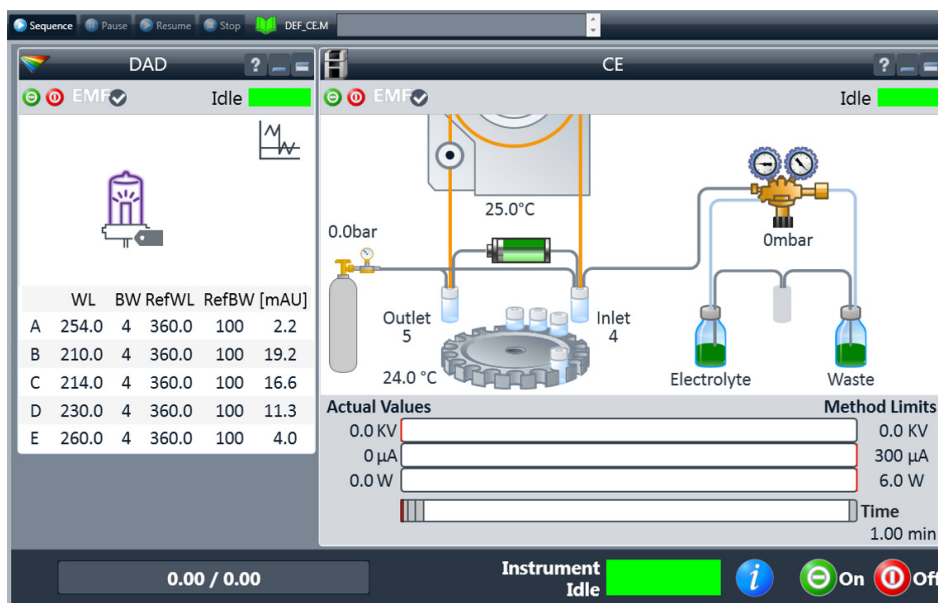
Figure 50 CE diagram with CEC mode activated

CEC-Specific GUI Explanation

When the **CEC mode** is selected, a new icon in the form of a gas cylinder appears on the left right side of the CE diagram next to the tray (see [Figure 50](#) on page 173).

There are also a number of new parameters which may be set either from the **Edit Entire Method** option or from the **Individual Parameter Settings** within the **Instrument** menu selected from the **Method & Run Control** toolbar. The new parameter choices appear in the following menu selections.

CEC Method Setup

When **CEC mode** is selected, a new section for pressure settings appears on the left side of the **Method Setup** dialog box. (see [Figure 51](#) on page 174). The parameter set here is the pressure applied to both vials in the inlet and outlet lift during the electrophoretic run.

Options that can be selected include:

5 Using the Agilent 7100 Capillary Electrophoresis System Capillary Electrochromatography, CEC

- Pressure off
- Internal, pressure or vacuum up to 100 mbar can be applied to the inlet lift
- Raw External Pressure
- External, set pressure from 2-12 bar

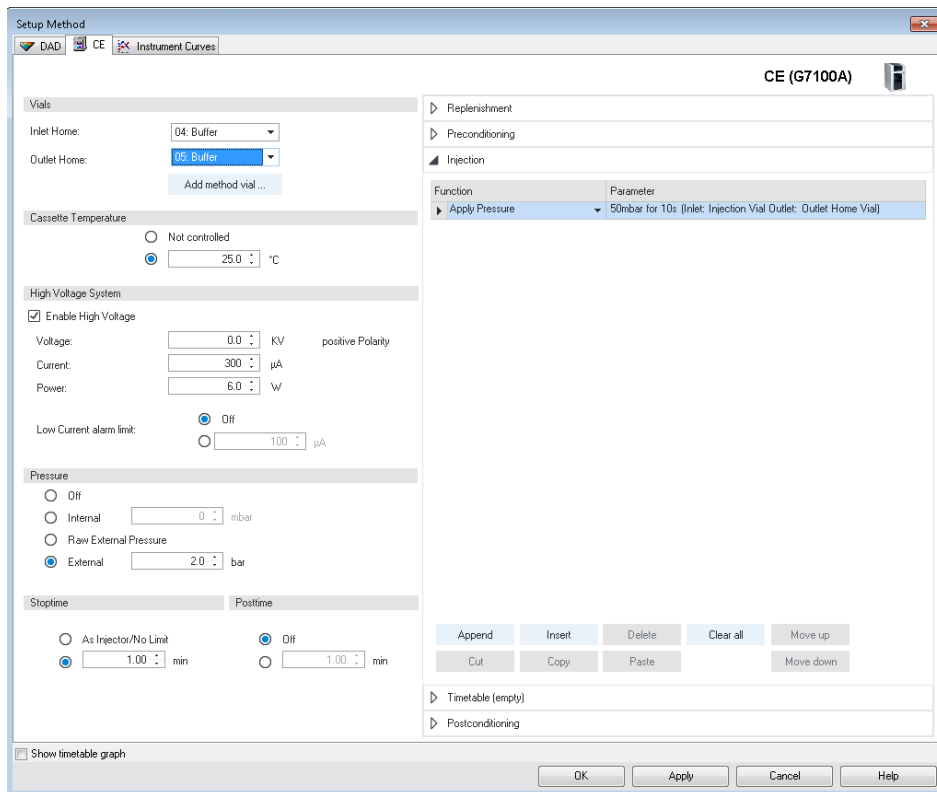
Figure 51 CEC setup method screen

Preconditioning and Postconditioning

In order to condition a packed capillary, an applied pressure of approximately 12 bar alone is usually insufficient to move the mobile phase rapidly through the column. Therefore, the applied high pressure may be assisted by the application of an electric field. This augments the high pressure flow with an electroosmotic flow. In the preconditioning or postconditioning section of the set-up method, the **High pressure flush** command can be used for this purpose, see [Figure 52](#) on page 175.

High pressure can be applied to the capillary inlet vial which may be:

- A set value (2-12 bar)
- Raw external pressure (up to 15 bar): Allows pressurizing the inlet vial and/or the outlet vial. The external pressure up to 15 bar can be applied directly without reduction.

High pressure flush

Raw Pressure
 bar

KV

μ A

W

s

Inlet

Outlet

Apply pressure to Outlet

Apply pressure to both

Figure 52 High pressure flush command menu

5 Using the Agilent 7100 Capillary Electrophoresis System

Capillary Electrochromatography, CEC

Other parameters which can be set in the **High pressure flush** command menu include:

- Applied voltage (kV), current (uA) or power (W)
- Duration (s)
- Inlet vial and/or outlet vial

NOTE

When applying voltage simultaneously, the inlet and outlet vials must contain electrolyte in order to maintain electrical contact.

Injection Options with CEC

The installed high pressure supply may also be used for injection purposes using the **Timetable** option in the **Injection** field of a method setup.

From this menu you may select the **High pressure flush** command during injection with the same parameter choices as those available during preconditioning.

This may be used to apply a pressure injection while using a replaceable gel-filled capillary.

NOTE

Please note that the external pressure is not as precise as the 50 mbar injection pressure!

Timetable Options with CEC

The installed high pressure supply may also be used in the general timetable of a method setup. The **Pressure** command is extended by the **External** pressure functionality, see [Figure 53](#) on page 177.

Figure 53 Pressure options in CEC mode

Timetable-controlled changes in the pressure settings are:

- OFF: Pressure usage is switched off.
- Internal, pressure (-100 mbar to +100 mbar): Can be applied to the inlet vial.
- Flush: Standard flush procedure is applied with an internally generated pressure of approximately 950 mbar.
- Raw external pressure (up to 15 bar): Allows pressurizing the inlet vial and/or the outlet vial. The external pressure up to 15 bar can be applied directly without reduction.
- External: Allows pressurizing the inlet vial and/or the outlet vial. External pressure is reduced to the set value between 2 and 12 bar and applied in a controlled manner.

CEC Control Using the CE Diagram

High pressure may be applied independently of the method for capillary conditioning or cleaning by clicking right at the external pressure icon in the CE diagram (see [Figure 50](#) on page 173). This will cause a parameter setting field to appear (see [Figure 51](#) on page 174). From this field you may chose to apply pressure to the inlet, outlet or both vials, and also whether to simultaneously apply a voltage of -30 to 30 kV for a given time.

Running the CEC Analysis

After setting up the operational parameters, you can initiate the run from the **Method & Run Control** menu by clicking **Run Method** or by clicking **Single Sample** in the system diagram. Some other operational aspects should be noted when running CEC separations.

- Capillaries** Packed capillaries are very fragile and must be handled with great care. Offline purging of the capillary using an LC pump may be necessary if air bubbles form within the capillary.
- Buffers** CEC buffers generally have lower concentrations (4–20 mM) than conventional CE buffers. These may also contain a larger proportion of organic modifier than used in CZE or MEKC buffers. The currents generated are generally on the order of 2–10 μA for an applied field of approximately 600 V/cm.
- Injection** Because of the resistance of the packed bed, the usual applied pressure of up to 50 mbar is insufficient to introduce sample onto the capillary. In this case, high pressure and/or electrokinetic loading may be used, although the limitations of this technique should be noted.

Problem Solving


- Offline purging** If air bubbles occur during an analysis, the current will become instable or very low ($< 0.3 \mu\text{A}$). If this occurs, remove the packed capillary and connect it to an LC pump capable of delivering a flow rate of 40 $\mu\text{L}/\text{min}$ to the capillary.

NOTE

Great care should be taken when removing and handling the capillary and when connecting and disconnecting the capillary to or from the pump. Fused silica packed capillaries are extremely fragile.

- Instrument purging** If the capillary is filled with solvent, then this may be replaced by inserting the capillary into the instrument and flushing with external pressure using the external gas supply icon on the CE diagram. Generally, 10–60 minutes are sufficient to replace the storage solvent with the operating mobile phase, depending on the velocity of EOF.

5 Using the Agilent 7100 Capillary Electrophoresis System
Capillary Electrochromatography, CEC

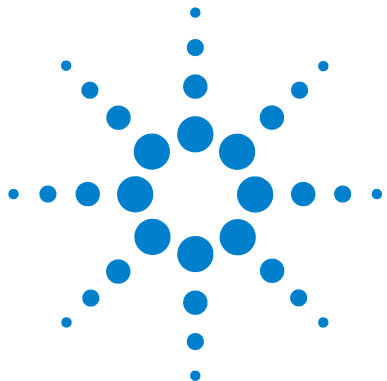6 Troubleshooting and Diagnostics

CE Problem Solving 182

Agilent Lab Advisor Software 189

This chapter gives an overview of troubleshooting and diagnostic features and the different user interfaces.

CE Problem Solving

The problem solving section in the instrument online help assists you in solving the most common failures. Nevertheless, a short list of problems is featured here.

Power On Fail

The instrument will boot up after turning the instrument on by pressing the power on switch. A green LED in the power switch is illuminated. If the instrument power on fails, the green control LED in the power switch may not light up.

- Check the external power supply, power cords and plugs for proper connection.
- Check the line fuses next to the power plug on the right rear of the instrument.
- Check fuses 1 to 6 on the main electronic board. A blown fuse is indicated by a small red LED visible through a hole in the metal shield of the instrument and a flashing green LED at the power switch.

Leaks

The leak sensor in the Agilent CE instrument is located in the lower left part of the instrument, see [Figure 54](#).

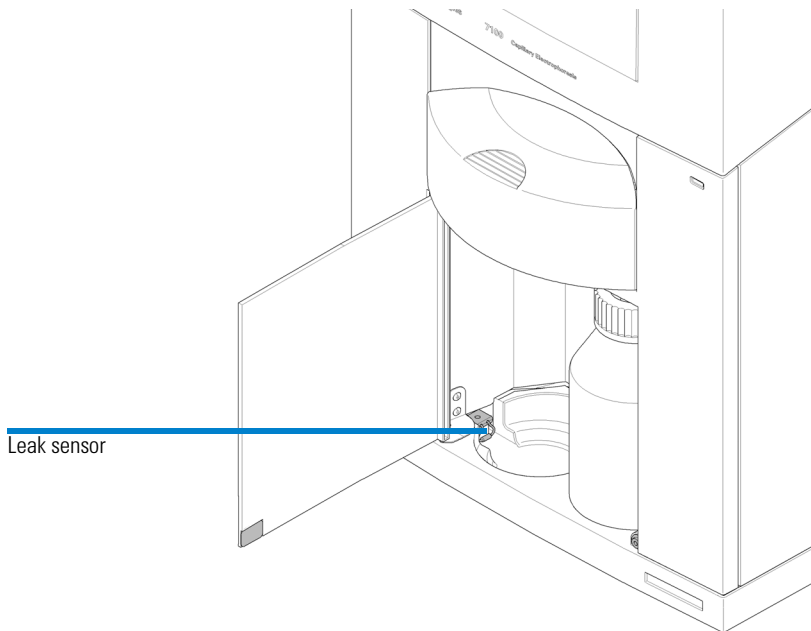
Figure 54 Leak sensor position

When the leak sensor is activated:

- The error LED on the front of the instrument is illuminated.
- The CE status displays an error.
- A logbook entry is created.
- The run is stopped.
- Pressure is released.
- High voltage application is stopped.

If the leak sensor is activated or you think there is a leak in the instrument:

- 1 Immediately switch off the Agilent CE instrument (and the external water bath, if connected).
- 2 Disconnect the line power cord.
- 3 Open the lid.
- 4 Remove the front cover as described to diagnose the cause of leak.
- 5 Dry the leak sensor using a tissue.

NOTE

The error will stay active for the ChemStation or Lab Advisor even after corrective action and drying. Please acknowledge the error by clicking "Clear error" in the Lab Advisor or click on the green poser button in the top left corner of the CE system diagram in the ChemStation.

Possible Reasons for Leaks and Appropriate Action for Spilled Liquid from Other Sources

- If you have spilled some liquid into the Agilent CE instrument, it is sufficient to thoroughly clean and dry the Agilent CE instrument and the leak sensor.

Leaking Water Bath Tubing Inside the Agilent CE Instrument

- Call Agilent Technologies.
- Do not use the Agilent CE instrument until Agilent Technologies has repaired it.

Leaks in Replenishment Tubing or Valves

- Call Agilent Technologies.

Reason Unknown

- If you can not trace the reason for the leak, call Agilent Technologies. There may be an instrument problem.

Blocked Replenishment Needle

If you are using the replenishment function of the Agilent CE instrument, buffer may precipitate inside the replenishment needle. To avoid this, put a vial filled with water or buffer (to a height of 1.8 cm) in vial position 49. Vial position 49 is the home position of the replenishment lift. The vial will be picked up by the lift and brought to the replenishment needle. The needle will dip into the water, thus preventing the buffer from precipitating. (see [“Preparing Vials for Buffer and Sample”](#) on page 129, Vial position 49)

Leakage Current

Leakage current is indicated by a yellow bar in the current part of the energy status display in the CE diagram. Only high leakage current will lead to this. An instrument curve allows you to monitor current leakage even at low values from the ChemStation.

Leakage current means that the current at the inlet electrode is different from the one at the outlet electrode and thus a current flows outside the capillary. A certain leakage current must be regarded as acceptable, since it is not avoidable due to technical reasons.

It can be caused by:

- A broken capillary
- An empty injection vial
- An air bubble in injection vial/or run buffer vials
- An air bubble in capillary
- A contaminated insulation plate
- A high humidity level
- A conductive capillary surface

If you find a broken capillary in the alignment interface, use an ultrasonic bath to clean the interface and remove quartz residues.

Broken Capillary

To check for a broken capillary, remove the capillary cassette and open it as described under “[Removing the Capillary Cassette from the Agilent CE Instrument](#)” on page 135. The capillary is especially sensitive to breaking at the detection window and the capillary ends, therefore check these parts of the capillary. If the capillary is broken at the detection window, a leakage current of nearly the same value as the expected current may occur.

Small instabilities in the current curves, as shown in [Figure 55](#), may also be a sign of a broken capillary.

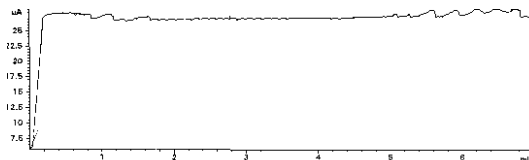
Figure 55 Current instabilities

If the capillary is broken at the ends, it may no longer reach into the liquid in the vial. In that case, the analysis is stopped if the low current alarm limit has been reached (this must be set first). Sometimes the capillary has very small cracks which are hardly recognizable under a microscope.

- Exchange the broken capillary

Contaminated Insulation Plate

Salt crystals or other contamination of the insulation plate may cause leakage current.

- Clean the insulation plate as described in “[Cleaning the Insulation Plate](#)” on page 227.

Empty Injection Vial

Access the vial position specified in the **Sample info** or the **Sequence table** using the **Get vial** function in the **Tray control** screen.

- Check if the vial is filled.

Dirt in the Vials

Clean all vials prior to use. First, flush them with water followed by, for example, some isopropanol.

- Use compressed air to dry the vials inside.

Air Bubble in Injection Vial or Run Buffer Vials

When filling sample or buffers into the vial, an air bubble is sometimes formed. This happens very frequently when using microvials. Air bubbles can cause air to be injected instead of the sample.

- To avoid air bubbles, carefully fill vials starting from the bottom.
- If you cannot avoid the formation of air bubbles, adjust the lift offset in the **Home values** dialog box so that the capillary doesn't reach into the bubbles.

Air Bubble in Capillary

Small air bubbles in the capillary may cause spikes in your electropherogram. Larger air bubbles expand so much during analysis that the current drops below the lower alarm limit.

- Flush your capillary with buffer to remove air bubbles.

Problems with the Replenishment System and Buffers Containing SDS

When a vial is replenished the SDS foam may stick to the side of the vial. The empty function draws air around the foam into the replenishment needle. When filling this vial, a set of new foam bubbles is introduced into the vial before the fresh buffer. The surface sensing system recognizes the first foam bubble as liquid. So the final liquid level is not at the height it should be. It may also happen that the vial contains no buffer liquid but only foam. The result could be seen as low current or even no current supplied.

As a workaround we recommend programming a replenishment program in two steps.

- 1 Replenish to a level about 0.2 cm higher than required.
- 2 Fill to the desired level.

If the fill results a level higher than the desired one, the vial is emptied to the level without using the sensing function and the foam has no influence.

As an example: the desired height is 1.4 cm

- 1 Replenish vial XX to 1.6 cm (0.2 cm more)
- 2 Fill vial XX to 1.4 cm

Alternative: No Replenishment Necessary.

- 1 Fill several vials with buffer.
- 2 Program methods which use these vials as the inhome and outhome vials (for example vials 4, 5, 6, 7).

Method 1:

inhome = 4, outhome = 5

Method 2:

inhome = 6, outhome = 7

Agilent Lab Advisor Software

The Agilent Lab Advisor software is a stand alone product that can be used independently of a data system. Agilent Lab Advisor software helps to manage the lab for high quality electrophoresis and chromatographic results and can monitor in real time a single Agilent CE or all the Agilent GCs and LCs configured on the lab intranet.


Agilent Lab Advisor software provides diagnostic capabilities for all Agilent 7100 CE instruments. This includes diagnostic capabilities, calibration procedures and maintenance routines for all the maintenance routines.

The Agilent Lab Advisor software also allows users to monitor the status of their CE instruments. The Early Maintenance Feedback (EMF) feature helps to carry out preventive maintenance. In addition, users can generate a status report for each individual CE instrument. The tests and diagnostic features provided by the Agilent Lab Advisor software may differ from the descriptions in this manual. For details, refer to the Agilent Lab Advisor software help files.

Configuration

The Lab Advisor software can communicate with a 7100 CE instrument via LAN or by a serial connection (RS232 cable, null-modem). To configure a connection, go to the section Configuration in the Lab Advisor software. Select **Instrument > Available Instruments** and as the instrument type **Agilent CE**. You can enter either the LAN address (for example 192.168.254.11) or the respective COM port of your PC (for example Com1) under instrument address.

6 Troubleshooting and Diagnostics
Agilent Lab Advisor Software

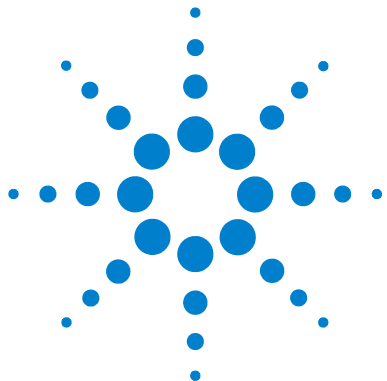7 Hardware Information

Identifying Instrument Components	192
7100 Capillary Electrophoresis Instrument Electronics	193
Firmware	196
Diode Array Detector	199
Changing the Fuses for the Power Supplies	202
Installing Drainage Tubing	203
External Water Bath for Tray Cooling (Optional)	204
Analog to Digital Converter	206

This chapter provides a more detailed description of the instrument's hardware and electronics.

Identifying Instrument Components

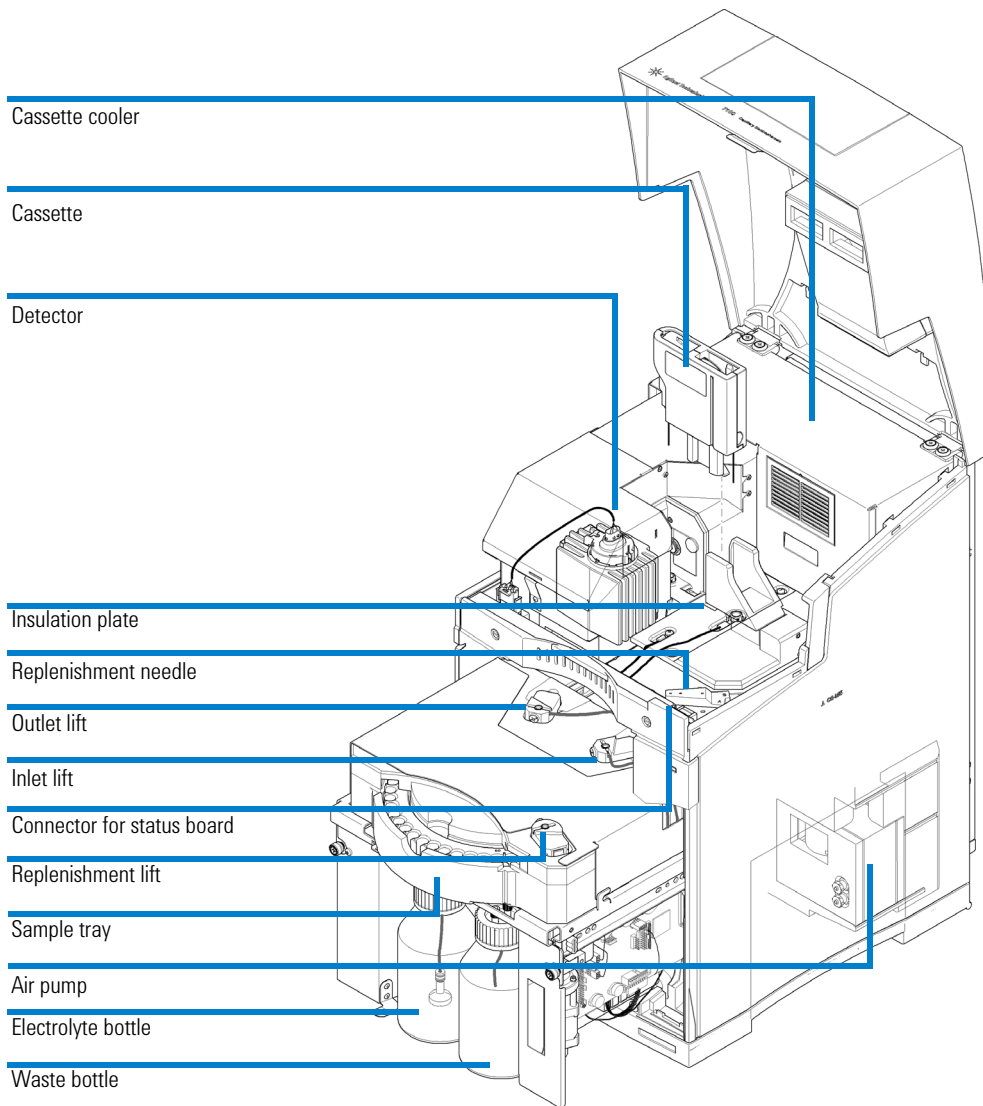
Figure 56 Overview of instrument components

7100 Capillary Electrophoresis Instrument Electronics

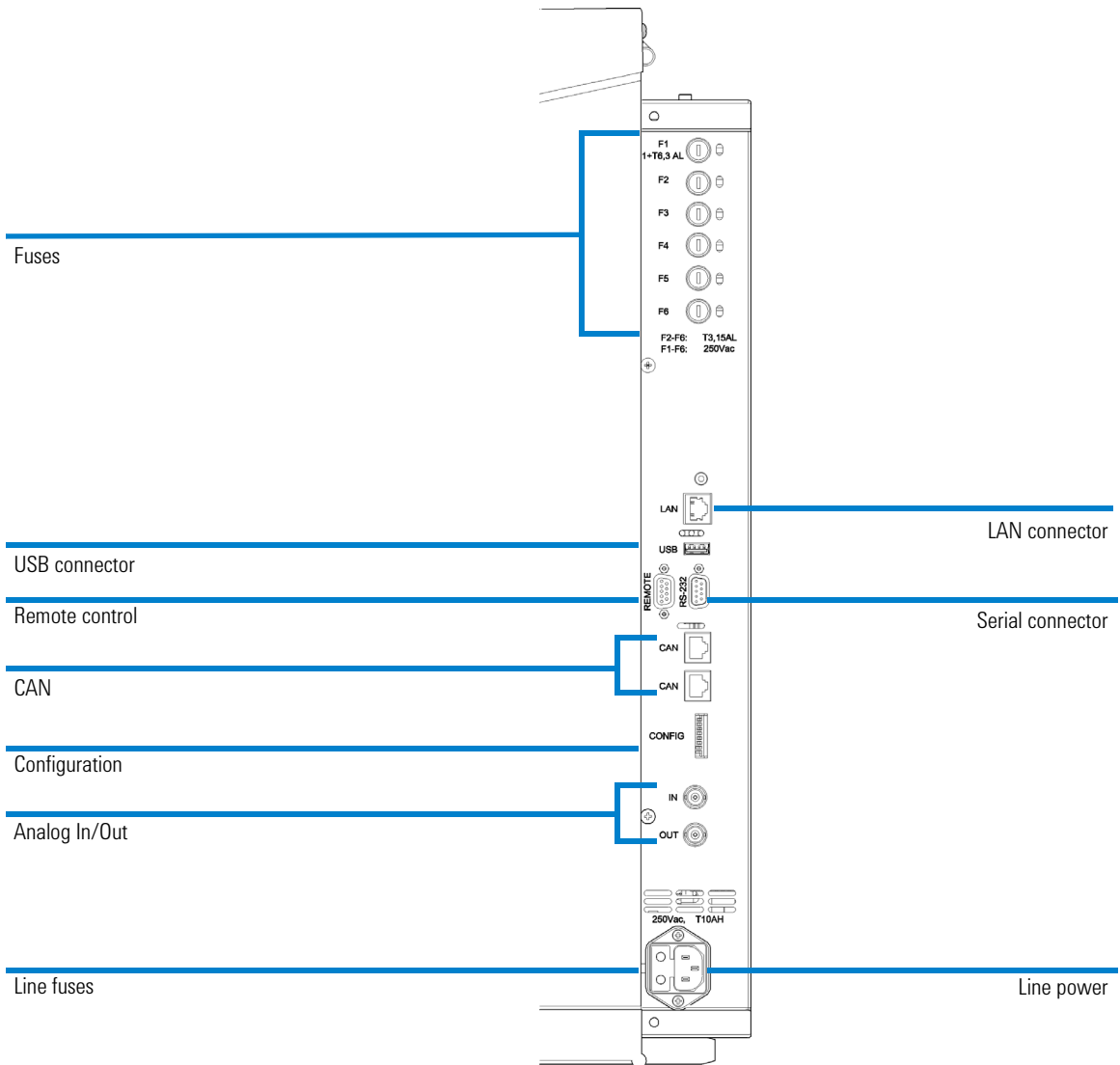
Figure 57 Connectors on the rear of the instrument

All electrical connectors are placed on the right side of the instrument, see [Figure 57](#) on page 193.

Available Connectors:

- Power connector, Fuse netfilter: 2 fuses, 250 VAC, T6.3AH, (**part number 2110-1018**)
- LAN connector (7100 to controlling PC)
- CAN connectors (7100 to additional Agilent 1200 Series Module)
- USB connector (for future use)
- RS232 connector (serial interface to a computer. This connection has a special out pin and is not compatible for connecting printers or plotters. It is called "null modem cable" with full handshaking and the wiring is established between pins 1-1, 2-3, 3-2,4-6, 5-5, 6-4, 7-8, 8-7 and 9-9.)
- APG remote connector (Agilent Technologies Analytical Products Group remote connector to be connected to Agilent modules, for example to MS system)
- Analog out (BNC connector, 0 to 0.1 V or 0 to 1 V)
- Analog in (BNC connector, 0 to 1 V, shield is analog - and center is analog +)
- 8-bit configuration switch (see [Chapter 4](#), "LAN Configuration," starting on page 75)

Fuses

- Fuse 1 (Peltier) T6.3AL, 250 VAC, (**part number 2110-0623**)
- Fuse 2 (HV supply, CE detector board) T3.15 AL, 250 VAC, (**part number 2110-1417**)
- Fuse 3 (CE sampler board with valve and lifter control) T3.15 AL, 250 VAC, (**part number 2110-1417**)
- Fuse 4 (heater, optional), T3.15 AL, 250 VAC, (**part number 2110-1417**)
- Fuse 5 (processor, +5 V, +15 V, -15 V supplies on mainboard, air pump, fans, CE temperature sensor board and cassette fan) T3.15 AL, 250 VAC, (**part number 2110-1417**)
- Fuse 6 (DAD) T3.15 AL, 250 VAC, (**part number 2110-1417**)

NOTE

Next to each fuse F1 to F6 is a LED on the rear of the electronic board. This is visible through a cutout in the metal shield of the electronic box. A red LED indicates that the respective fuse is blown. If one of the fuses is blown, the green LED of the power switch flashes.

WARNING**Fire hazard**

Using wrong fuses may result in a fire hazard.

- For continued protection against fire hazards, only replace fuses with the same type and rating.
 - Only use the fuses specified for this instrument. Use of other fuses or materials is prohibited.
-

Firmware

The firmware for the instrument consists of two independent sections:

- A non-instrument-specific section, called the resident system.
- An instrument-specific section, called the main system.

Resident System

The resident section of the firmware is identical in many Agilent devices and modules. Its properties include:

- The complete communication capabilities (CAN, USB, LAN and RS-232C)
- Memory management
- Ability to update the firmware of the main system

Main System

Its properties include:

- The complete communication capabilities (CAN, USB, LAN and RS-232C)
- Memory management
- Ability to update the firmware of the resident system

In addition, the main system comprises the instrument functions that are divided into:

- Common functions:
 - Run synchronization through APG remote
 - Error handling,
 - Diagnostic functions
- Or module-specific functions:
 - Internal events such as lamp control, filter movements
 - Raw data collection and conversion to absorbance

Firmware Updates

Firmware updates can be done using the Lab Advisor Software, PC and Firmware Update Tool with local files on the hard disk.

To replace or update the firmware:

- 1 Connect to the instrument with Lab Advisor and refer to the section **Firmware Update**.
- 2 Provide the firmware file on your local hard disk. You might locate the file in any folder and provide the location by selecting **browse to folder** or put it to a folder specific for your instrument (CE/LC).
- 3 Connect to the instrument, select the file in a standard folder or browse to the respective folder. Download it to the instrument.

NOTE

While the firmware is downloaded to the instrument, the status indicator LED on the front of the instrument front will blink in red.

- 4 Disconnect the firmware update tool from the instrument.

You can find a detailed description in respective help files of the Lab Advisor.

The file naming conventions are:

PPPP_RVVV_XX.dlb (for example **7101A_U.06.23_0002.dlb**).

PPPP represents the product number, R the firmware revision, VVV is the revision number and XX is the build number of the firmware.

NOTE

Updates to main system can only be done in the resident system. Updates to the resident system can only be done in the main system.

Main and resident firmware must be from the same set (please refer to the Agilent HPLC Firmware Bulletin for more details).

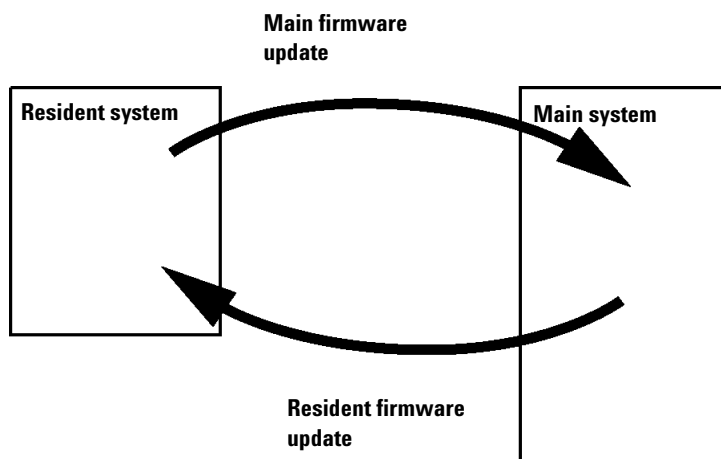
Figure 58 Firmware update mechanism

The latest firmware revisions and documentation are available on the Agilent website (<http://www.chem.agilent.com>)

Diode Array Detector

The DAD is directly controlled by the DAD diagram, through method parameters and general control settings.

Direct control of the DAD and control through method parameter is described in Controlling the Agilent CE instrument (“[Detector Parameters](#)” on page 121) and the section “User interface” in the ChemStation online help.

DAD Control

In the **Method and Run Control** of the ChemStation select **Instrument > More DAD > Control... > Control** to open the DAD control dialog (Figure 59 on page 200).

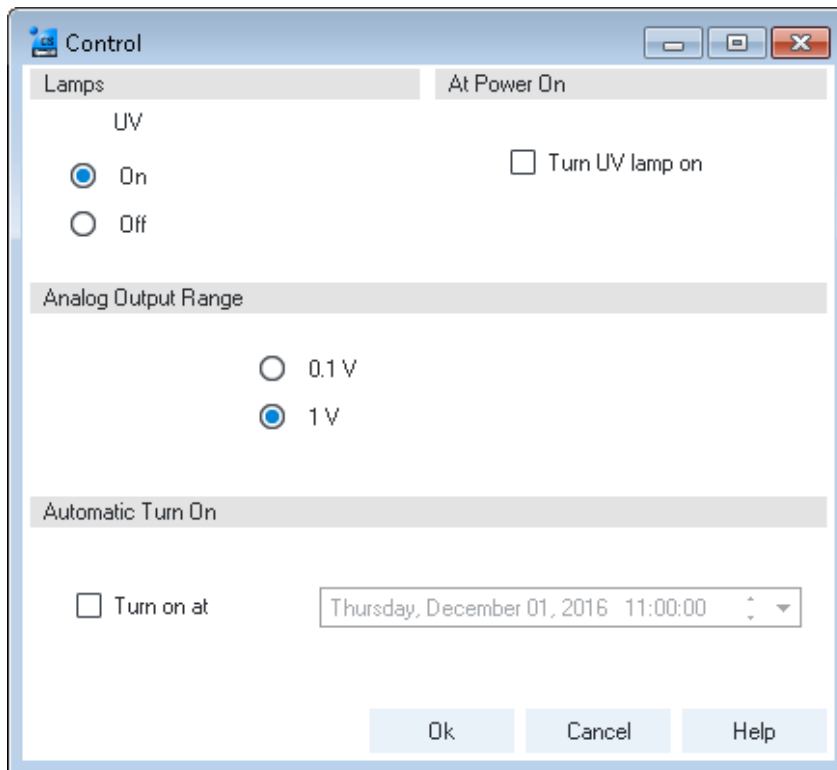
Figure 59 DAD control dialog

You may switch the UV lamp on or off in the **Lamps** dialog. Changes are applied only after clicking **OK**. A checkmark in the section **At Power On** allows you to immediately turn on the deuterium lamp when the instrument is switched on.

In the **Automatic Turn On** field you can also set a date and time when the lamps are switched on automatically. The instrument must be kept powered on, for example overnight, to use this feature.

Digital to Analog Converter

The 7100 CE System provides a direct analog output of chromatographic signals A. It may be used to connect to another data acquisition system or chart recorders. The analog signal is converted from digital format by the on-board 20-bit digital-analog-converter (DAC). Please select the suitable voltage range for your analog recording hardware under **Analog Output Range**.

Changing the Fuses for the Power Supplies

Prior to changing fuses, you must switch the instrument off and disconnect it from the line power.

Fuses are located on the back right side of the instrument. Replace fuses with identical part numbers as specified under “[7100 Capillary Electrophoresis Instrument Electronics](#)” on page 193.

WARNING

Fire hazard

Using wrong fuses can result in fire hazard.

- For continued protection against fire hazard, replace fuses only with the same type and ratings.
 - Only use the fuses specified for this instrument. The use of other fuses or materials is prohibited.
-

Installing Drainage Tubing

The tubing drains condensation out of the instrument which may appear if you connect an external water bath to cool your samples in the sample tray.

On the left rear side of your instrument you may install silicone tubing. Place a beaker of at least 200 mL capacity under this tubing to collect the drained condensation.

External Water Bath for Tray Cooling (Optional)

The two connections on the left rear of the instrument (see [Figure 2](#) on page 73) can be used to install an external water bath for the tray of the Agilent CE instrument.

The temperature for the cooling liquid is set at the external water bath. The temperature display in the CE diagram displays the actual temperature of the tray controlled by the cooling liquid. The external water bath cannot be controlled by the Agilent CE instrument.

Good Practices

To avoid problems when using the external water bath, consider the following points.

- The tubes used should have thermal insulation to avoid condensation.
- Use a displacer to minimize the amount of liquid that is released in case of leaking cooling lines.
- The device should be equipped with a liquid level sensor to switch off automatically if liquid is lost.
- Do not operate the devices at temperatures below +1 °C to avoid formation of ice on internal surfaces.

External Water Bath Requirements

- Minimum cooling power 100 W, at bath temperatures of 5 °C.
- Temperature precision better than ± 0.2 °C.
- Safety requirements must meet national standards and also: be suitable for unattended operation, be suitable for continuous operation, and have controllable high temperature protection.
- Requirements for external liquid circulation:
 - Suitable for external circulating of liquid,
 - Suitable for connection of 8 mm ID tubes,
 - Minimum circulating liquid flow rate 8 L/min,
 - Minimum differential pressure of 150 mbar, and
 - Maximum pressure of 500 mbar.

- Requirements for pressure pump:
A pump with the capacity to generate a differential pressure of at least 0.15 bar. For a single action pump, a high pressure capability of 0.15 bar is required. The high pressure must not exceed 0.5 bar and the low pressure must not exceed -0.5 bar.
- Cooling liquid:
Distilled water plus a maximum 3 % anti-freeze.
- Bath temperature:
The minimum allowed temperature is 1 °C and the
Maximum allowed temperature is 50 °C.

NOTE

While setting up the bath temperature, monitor the tray temperature in the user interface of the data system and make sure that the tray temperature is within the specification (+10 °C - +40 °C, see performance specifications) to avoid condensation and potential damages to the instrument.

Connecting an External Water Bath

- The startup kit supplied with the Agilent CE instrument includes two fittings for flexible tubing with 8 mm internal diameter. Use those fittings for the flexible tubing used with your external water bath.
- Connect the fittings to the connectors on the instrument (see [Figure 2](#) on page 73). Use the upper connector for incoming liquid, the lower connector for outgoing liquid.

NOTE

Make sure both the water bath and the Agilent CE instrument are switched off when installing the external water bath.

CAUTION

Risk of liquid spill and instrument damage

Higher pressure might lead to liquid leakage inside and outside the instrument. Elevated temperatures might lead to deformation and leakage of the tubing system. Liquid spillage can damage electronic components of the instrument.

- Do not use a water bath with more than 0.5 bar pressure.
- Do not use liquids at temperatures above 50 °C.

Analog to Digital Converter

Some detectors (for example CCD detectors) require an intermediate interface to convert analog data into a digital format that can be sent to the ChemStation (Analog In). The Agilent 7100 Capillary Electrophoresis System has a 20-bit AD converter built in. It enables voltage conversions ranging from 0 V to 1 V.

You can set up name, units and conversion factors for the analog in signal from the following dialog which is accessible under Instrument/configuration, CE-part configuration, see [Figure 60](#).


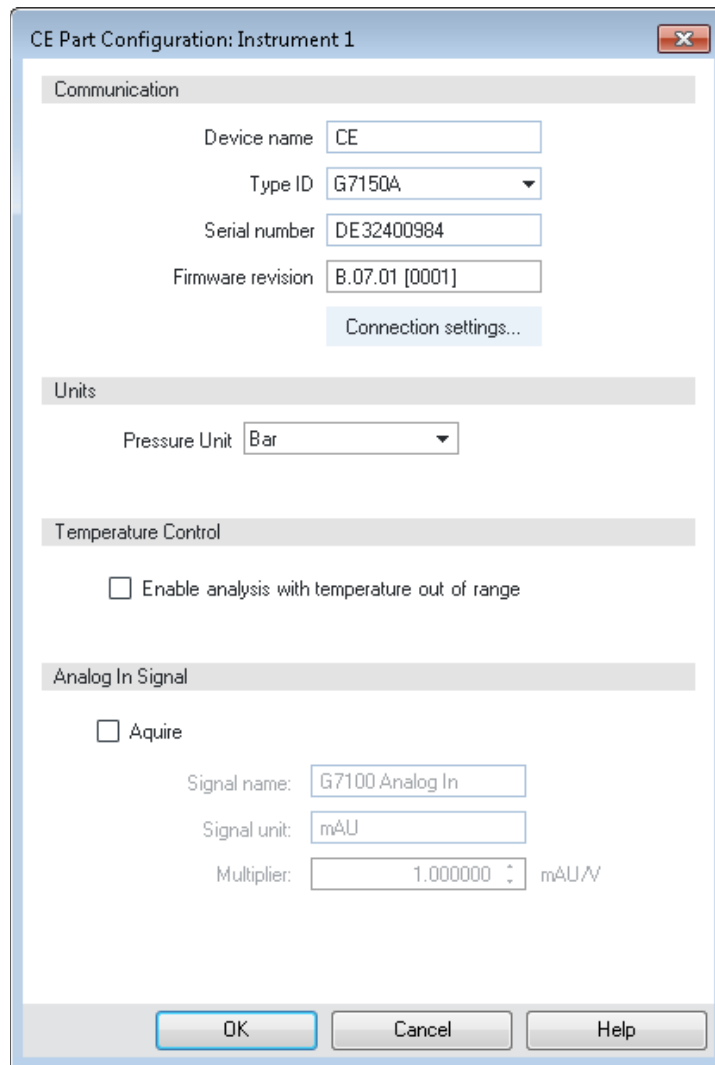
Figure 60 CE part configuration

7 **Hardware Information**
Analog to Digital Converter

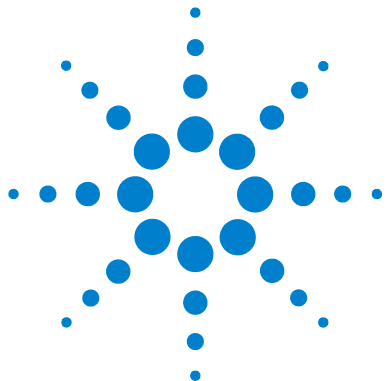8 Maintenance

Standard Operating Procedures for Maintenance of Your Agilent CE Instrument	210
Overview of Maintenance	211
Early Maintenance Feedback (EMF)	214
Cleaning the Electrodes, Pre-punchers and Insulation Plate	216
Maintenance of the Replenishment System	236
Changing the Air Inlet Filter	242
Changing the Lamp	244
Cleaning the Instrument	248

This chapter describes how to maintain the CE instrument.

Standard Operating Procedures for Maintenance of Your Agilent CE Instrument

This chapter describes standard operating procedures (SOPs) for regular maintenance of your instrument. Only perform maintenance on the Agilent CE instrument that is specified in this chapter. Other maintenance or repairs must be done by personnel trained by Agilent Technologies. Unauthorized maintenance may be dangerous and damages are not covered by warranty.

Maintenance tasks are supported by the Agilent Lab Advisor software or the Agilent Instrument Utilities software. It is delivered with the 7100 Agilent Lab Advisor software comes with a complete set of features. The Agilent Instrument Utilities software provides set of basic features.

These software packages assist you by providing hardware test suits, a set of tools used during procedures, Early Maintenance Feedback (EMF) parameters and scheduling functions. You can find a detailed description in the respective online help documentation.

Overview of Maintenance

Table 16 Overview of maintenance

Procedure	Typical frequency	Notes
Cleaning the electrodes, pre-punchers and insulation plate	Clean frequently*	<ul style="list-style-type: none"> Inspect pre-punchers for salt deposits, bent tips, deposits in top funnel. Replace if either electrodes or pre-punchers are bent, show signs of wear such as scratches, or cannot be cleaned.
Cleaning the electrodes	Clean frequently*	Inspect for salt deposits, O-ring integrity, bending, and check alignment.
Cleaning the detection window and the alignment interface		<p>Capillary window:</p> <ul style="list-style-type: none"> Gently clean with isopropanol and a lint-free cloth. <p>Optical alignment interface:</p> <ul style="list-style-type: none"> Inspect for occlusion under microscope or magnifier (dust or shards of polyimide). If capillary was broken while in the interface, inspect and remove fragments (ultrasonic device and air burst)
Preparing/cleaning the replenishment system	System cleaning should be performed when new buffer is added to the electrolyte reservoir.	Clean for storage with isopropanol/water. Filter all solutions prior to use (0.2 or 0.45 µm).
Maintaining the functionality		
Changing the air inlet filter		<ul style="list-style-type: none"> if the environment is very dusty if the system cannot generate enough pressure at least every three months.
Changing the lamp	Perform lamp intensity test from the detector test suite in the Lab Advisor regularly (for example bi-weekly) and monitor lamp life using EMF values.	Use the green-labeled alignment interface without capillary for tests.

8 Maintenance

Overview of Maintenance

Table 16 Overview of maintenance

Procedure	Typical frequency	Notes
Repair of replenishment needle		<ul style="list-style-type: none"> Use the Clean Level Sensor task to remove salt deposits inside the tubing system. Keep a water-filled parking vial in carousel position 49. Occasionally replace the parking vial in position 49 with water or buffer. Inspect for bending if clogged. Remove persistent clogging by a syringe flush with water/isopropanol or use ultrasonic device.
Buffer reservoir		<ul style="list-style-type: none"> Inspect O-ring in cap for damage, remove salt deposits. Do not use buffer reservoir for long-term storage (prevents oxygen saturation and bacteria from growing). Remove liquid from waste bottle.
Inlet frit		<ul style="list-style-type: none"> Inspect and clean (especially if absorbing buffers have been used). Replace when dirty.
Vial caps		<ul style="list-style-type: none"> Do not overuse vial caps (especially the PEO caps). Inspect for damage.

* The chosen frequency may be time-based, for example weekly or injection-based, for example 100 injections. It depends on your individual application and needs. The time and injection-based EMF counter can be customized.

NOTE

Do not open the top cover or remove the cartridge without lowering the lifts. Use the function maintenance position from the tools section of the Lab Advisor. Careless opening may lead to floating of electrodes with buffer or sample which leads to contamination or carry-over.

List of Standard Maintenance Procedures:

- Clean the electrodes
- Clean the replenishment needle
- Clean the pre-punchers
- Check alignment of electrodes
- Clean the replenishment tubing system
- Change the air inlet filter
- Change the buffer frit
- Change the UV lamp
- Clean the insulation plate
- Perform a (full) system diagnostic test

Early Maintenance Feedback (EMF)

Maintenance requires the exchange of components which are subject to wear or stress. Ideally, the frequency at which components are exchanged or cleaned should be based on the intensity of usage of the instrument and the analytical conditions, and not on a predefined time interval. The Early Maintenance Feedback (EMF) feature monitors the usage of specific components in the instrument, and provides feedback when the user-selectable limits have been exceeded. The visual feedback in the user interface provides an indication that maintenance procedures should be scheduled.

EMF Counters

The CE instrument provides EMF counters for the CE part and the detector. The counters increment with instrument use, and can be assigned a maximum limit which provides visual feedback in the user interface when the limit is exceeded. The counters can be reset to zero after the related parts are exchanged. The CE instrument provides the following EMF counters:

- Analysis start counter
- Accumulated air pump on time
- Accumulated UV lamp on time
- Number of ignitions

Using the EMF Counters

The user-settable EMF limits for the EMF counters enable the Early Maintenance Feedback to be adapted to specific user requirements. The useful lamp burn time is dependent on the requirements for the analysis (high or low sensitivity analysis, wavelength, etc.), therefore, the definition of the maximum limits need to be determined based on the specific operating conditions of the instrument.

Setting the EMF Limits

The setting for the EMF limits must be optimized over one or two maintenance cycles. Initially, no EMF limit should be set. When instrument performance indicates maintenance is necessary, take note of the values displayed by the instrument counters. Enter these values (or values slightly less than the displayed values) as EMF limits and then reset the EMF counters to zero. The next time the EMF counters exceed the new EMF limits, the EMF flag will be displayed, providing a reminder that maintenance needs to be scheduled.

NOTE

Please introduce self-made EMF counters to remind you to frequently clean the electrodes, pre-punchers and insulation plate.

Cleaning the Electrodes, Pre-punchers and Insulation Plate

Buffer precipitates, dust, moisture or other solid material in the electrodes and pre-punchers or under the insulation plate may cause arcing. They can lead to current leakage, buffer contamination or carry-over. Cleaning these parts regularly will prevent this.

The electrodes, pre-punchers and the insulation plate should be cleaned regularly. Depending on the application, cleaning frequency can be for example once a week, or if peaks show unusual tailing or unusual peak shoulders. Please be aware that the EMF counter analysis start lets you set customized limits for warnings. Cleaning is necessary if arcing or excessive current leakage is observed. Usage of creeping buffers, such as urea, may lead to higher cleaning frequency.

Steps for the cleaning procedure for electrodes, pre-punchers and the insulation plate:

- Prepare the Agilent CE instrument for this task
- Remove the covers
- Access the electrodes
- Access the liquid handling module
- Access the pre-punchers
- Clean the electrodes
- Clean the pre-punchers
- Clean the insulation plate
- Reinstall the pre-punchers
- Reinstall the insulation plate
- Reinstall the electrodes

NOTE

Follow the common safety regulations for laboratories when doing this task. Wear safety glasses and rubber gloves.

NOTE

Make sure you have disconnected the line power cord of the Agilent CE instrument before you proceed.

When	If arcing or excessive current leakage is observed.
Parts required	12 mm hexagonal socket screwdriver (included in the startup kit) Screwdriver for slotted screws (included in the startup kit) Screwdriver for cross head screws (included in the startup kit) Wash bottle with isopropanol Wash bottle with water Compressed inert gas, oil-free

Prepare the Agilent CE Instrument

Prior to maintenance do the following:

- 1 Make sure no analysis is running or will be started from the ChemStation.
- 2 Start monitoring your CE instrument with the Lab Advisor.
- 3 Select the **Tools** section in the instrument panel of the Lab Advisor.
- 4 Select the **CE control** screen and then **Maintenance positions**.
These steps unload all lifters and move them to the lowest position. With all lifters in this position, the tray is freely movable. A dialog reports which lifters were unloaded in case vials were in the system.
- 5 Turn off the Agilent CE instrument.
- 6 Disconnect the line power cord from the Agilent CE instrument.
- 7 Disconnect any of the following connection if present: waterbath, external pressure supply and drainage.

WARNING**Hot lamp**

If you were using the Agilent CE instrument shortly before, the lamp may be very hot.

- Avoid touching the lamp.
- Wear protective gloves.

Remove the Covers

- 1 Open the top cover.
- 2 Remove the capillary cassette (see “[Removing the Capillary Cassette from the Agilent CE Instrument](#)” on page 135)
- 3 Turn the front cover screws by a quarter of a turn to vertical orientation. They are not meant to be removed completely.
- 4 Open the door to the cabinet a little.
- 5 Carefully pull the front cover away from the main frame.
Pull steadily on both sides. You will need slightly more force to remove the right side due to a connector. Turn the screws back to horizontal orientation.

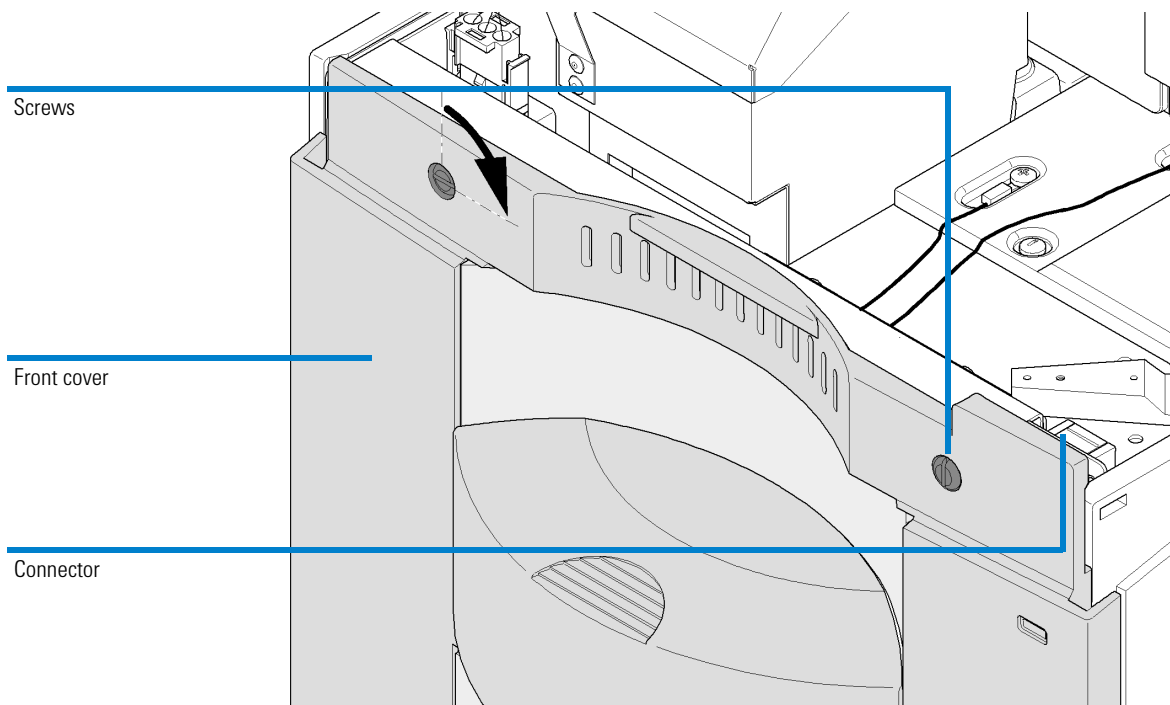
Figure 61 Removing the front cover

Accessing the Electrodes

- 1 Use a screwdriver to unscrew the two metal screws holding the insulation plate.
- 2 Use a screwdriver to remove the screw that secures the blue ground cable.

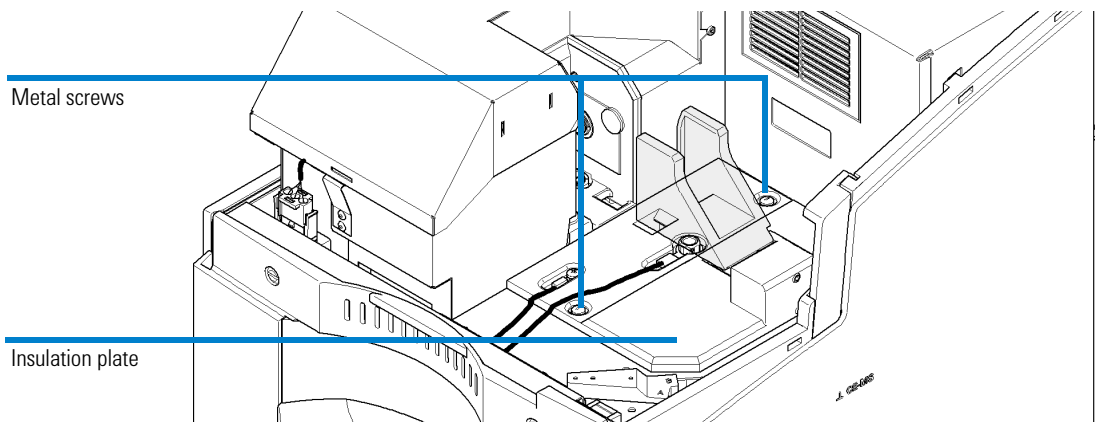
Figure 62 Insulation plate

- 3 Use the 12 mm hexagonal socket screwdriver to unscrew the inlet electrode.

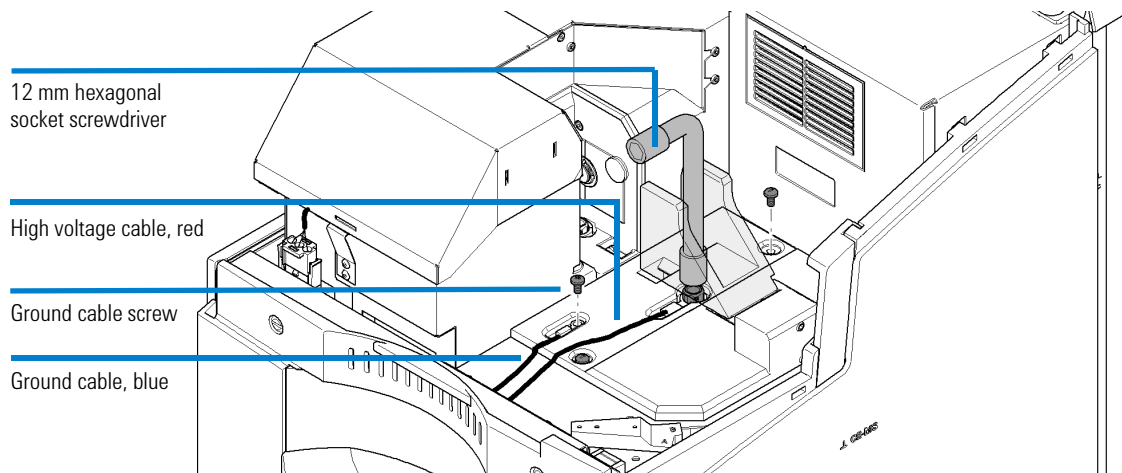
Figure 63 Removing the ground cable and the inlet electrode cable

8 Maintenance

Cleaning the Electrodes, Pre-punchers and Insulation Plate

- 4 Carefully lift the red high voltage wire. The inlet electrode will come out.
- 5 Slowly slide the insulation plate first to the right (1.), then tilt it up (2. and 3.) and finally take it out.

Make sure that the outlet electrode is not touching the foam of the tray cooling or the lift station.

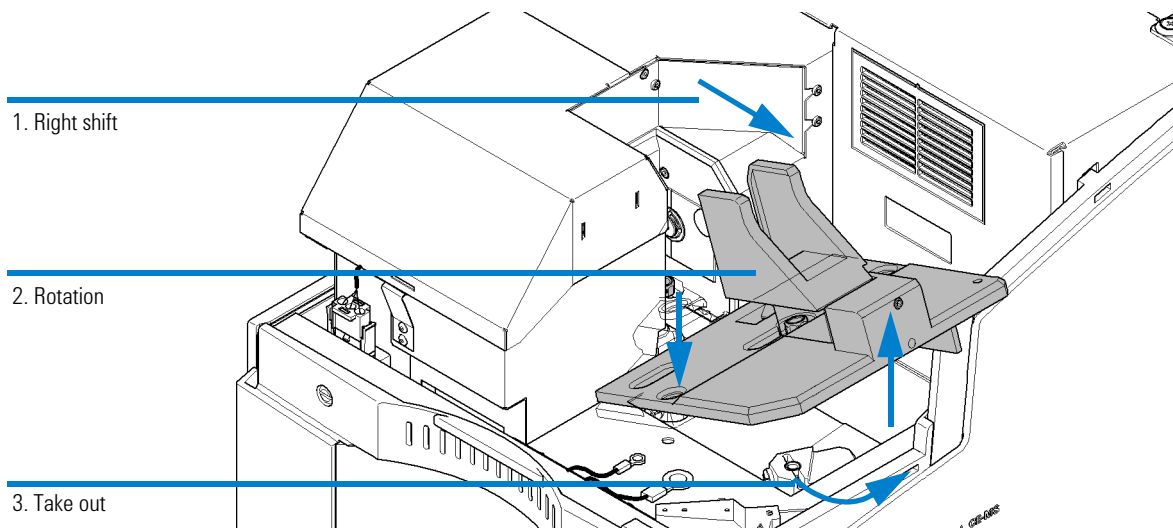
Figure 64 Removing the insulation plate

NOTE

It is not possible to remove the outlet electrode without taking out the insulation plate.

- 6 Remove the outlet electrode from the insulation plate using the 12 mm hexagonal socket screwdriver.

Accessing the Liquid Handling Module

CAUTION

Damage to tubes and liquid spills

Keeping tubes connected on the left side of the instrument while pulling out the liquid handling module will damage the tubes and might lead to liquid spills.

Disconnect any external pressure supply, drainage or connections to an external water bath for tray cooling prior to accessing the liquid handling module.

CAUTION

Damage to the instrument

Damage may lead to a non-operational instrument.

- Do not lift the liquid handling module upwards, outside the instrument. The module is connected to the main frame by two flat ribbon cables and two flexible tubes. When lifting it up, the liquid handling module might come off the guiding rail and the connections might be affected.
-

- 1 Use a screwdriver to unscrew the slotted screws holding the liquid handling module in the instrument.
- 2 Take the electrolyte and waste bottles out of the cabinet. Keep them connected and place them in front of the instrument.
- 3 Carefully pull the liquid handling module out from the main frame. Pull steadily in the middle until you reach the end position. Carefully avoid damaging the replenishment needle.

8 Maintenance

Cleaning the Electrodes, Pre-punchers and Insulation Plate

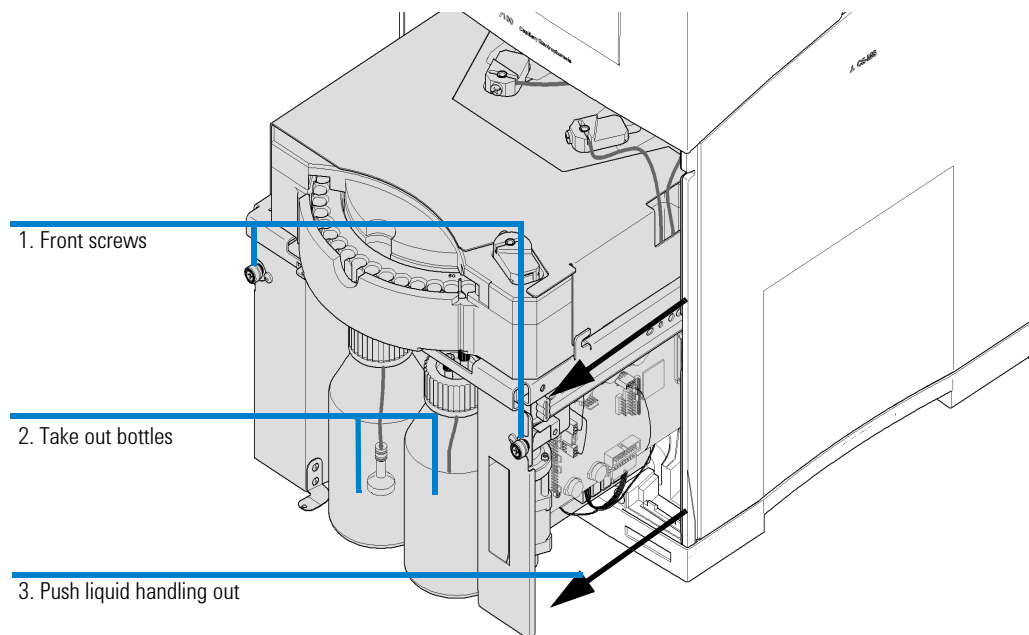
Figure 65 Accessing the liquid handling module

Accessing and Removing the Pre-punchers

Inlet and outlet lifts are located behind the sample tray in the rear of the liquid handling module. The replenishment lift is located in the front on the right side under the replenishment needle.

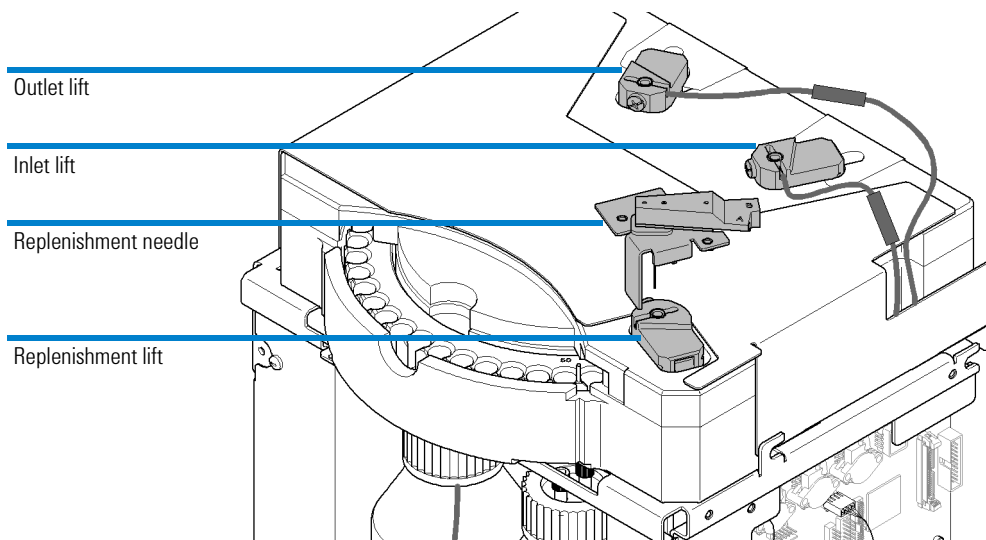
Figure 66 Inlet and outlet lifts, replenishment lift and replenishment needle

NOTE

Do not touch or bend the replenishment needle when handling the pre-punchers.

WARNING

Risk of injury due to sharp tips

The pre-punchers have very sharp tips.

- Be careful when touching the tips of the pre-punchers.

8 Maintenance

Cleaning the Electrodes, Pre-punchers and Insulation Plate

- 1 Remove the plastic screw which secures the pre-punchers using a Pozidriv screwdriver.

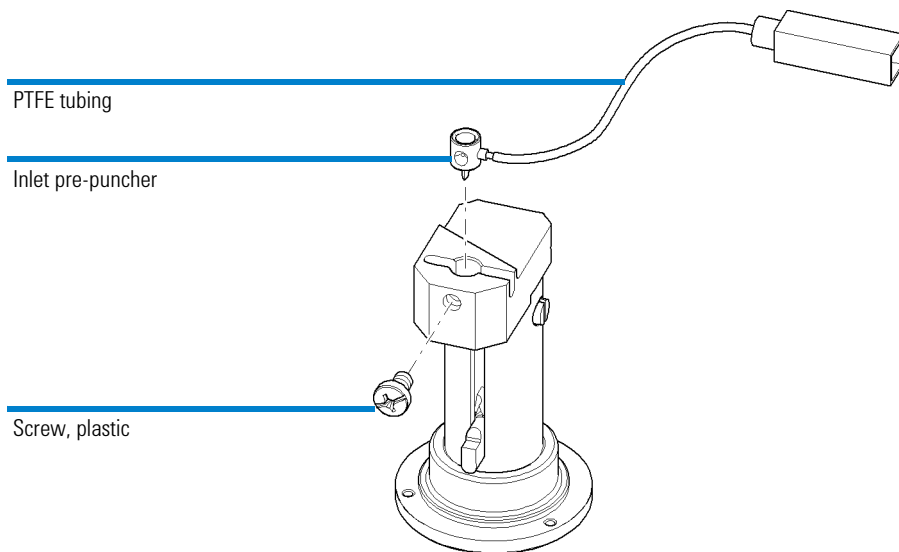
Figure 67 Plastic screw of the pre-puncher

- 2 The pre-puncher has a hole which can be used to gently lever it out using a paper clip. Carefully pull the pre-puncher out of the inlet, outlet and replenishment lift.
- 3 Unscrew the PTFE inlet pre-puncher tubing (labeled 7 on the tubing), the outlet pre-puncher tubing (labeled 8). The pre-puncher for the replenishment lift is not connected to a tube.

NOTE

The outlet lifter tubing (labeled 8) might be physically disconnected from the pressure system. This status is advantageous in case no high pressure applications are used. This connection is required for applications in CE+p and CEC mode.

NOTE

The PTFE tube does NOT need to be removed from the pre-punchers for cleaning. If the connection between the tube and the pre-puncher is not tight, since the tube became wider, shorten the tube and shift it on the pre-puncher again.

Cleaning the Electrodes

- 1 Flush the electrodes with water. Watch for salt crystals. All salt crystals must be removed.

Figure 68 Flushing electrodes

- 2 Flush the electrodes with isopropanol.
- 3 Use compressed air to dry the electrodes and remove all traces of liquid.

CAUTION

Increased leakage current and arcing due to wet electrodes

Increased leakage current and arcing might appear during analysis if electrodes are not completely dry.

- To avoid contamination, the inside of the electrodes must be completely dry before reinstallation.

NOTE

If there is still an accumulation of dirt in the upper funnel or inside the electrode start the cleaning procedure again. You may place it in a beaker of water or isopropanol and treat this in an ultrasonic bath for 5 minutes followed by steps 2 and 3 in this section. If this fails, replace the electrode.

Cleaning the Pre-punchers

- Check for damaged or bent tips before cleaning the pre-punchers.
If the tips are damaged, replace the pre-punchers.
- Check for grooves in the funnels of the pre-punchers.
Grooves are a sign of improper alignment of the electrodes and may lead to improper sealing. Improper sealing may lead to problems during flushing, injection or CEC applications. Use the alignment center of the Lab Advisor to align and check the installation. If the funnel shows grooves, replace the pre-punchers.

WARNING

Risk of injury due to sharp tips

The pre-punchers have very sharp tips.

- Be careful when touching the tips of the pre-punchers.

- 1 Flush the pre-punchers with water. Watch for salt crystals. All salt crystals must be removed. Take particular care with the PTFE tube connector.

Figure 69 Flushing pre-punchers

- 2 Flush the pre-punchers with water and isopropanol.
- 3 Use compressed air to dry the pre-punchers and remove all traces of liquid. Do not forget to dry the connected tubes as well.

NOTE

If there is an accumulation of dirt in the pre-puncher funnel, put the pre-punchers and connecting tubing in a beaker of water and treat it in an ultrasonic bath for five minutes followed by steps 2 and 3 in this section.

NOTE

Residual liquid in connecting tubes may be transported to the air valve blocks and result in damage. Such liquid plugs may also negatively influence injection reproducibility.

Cleaning the Insulation Plate

Salt crystals, films of substances such as buffer or gel that allow moisture will lead to increased leakage current or even arcing. The insulation plate must be free of such contamination.

- 1 Moisten a tissue with water and use it to clean the insulation plate.
- 2 Moisten a tissue with isopropanol and use it to clean the insulation plate.
- 3 Dry the insulation plate with a tissue and compressed air.

Reinstalling the Pre-punchers

- 1 Carefully insert the pre-puncher (labeled 8) in the outlet lift. Connect the tube to its connector (labeled 8) if you run applications in the CE+p and CEC mode. The fitting of the outlet pre-puncher must be tight. Otherwise, problems may occur during CEC mode.

NOTE

If you do not run applications in CE+p and CEC mode, keep the tube disconnected and fix the tubing in such way that it does not disturb the movement of the inlet lifter.

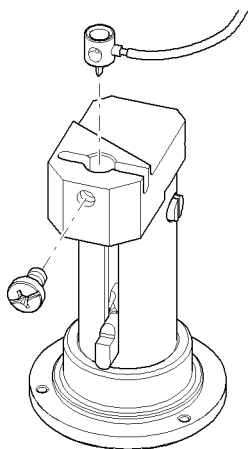
Figure 70 Inserting the pre-puncher

- 2 Carefully insert and tighten the plastic screw to secure the pre-puncher. The pre-puncher will assume the correct position as the screw is tightened.
- 3 Carefully insert the inlet pre-puncher with the air pressure tube (labeled 7) connected in the inlet lift. Connect the tube to its connector (labeled 7). The fitting of the inlet pre-puncher must be tight. Otherwise, pressure or injection problems may result.
- 4 Carefully insert and tighten the plastic screw to secure the pre-puncher. The pre-puncher will assume the correct position as the screw is tightened.

NOTE

The screws used for fixation of the inlet and outlet pre-puncher are made of plastic. Do not use metal screws. Metal screws may lead to leakage current and arcing which disturbs applications and may damage the instrument.

- 5 Carefully insert the replenishment pre-puncher. Tighten the screw to secure the pre-puncher.

The pre-puncher will assume the correct position as the screw is tightened. The replenishment pre-puncher is not connected to a tube.

Reinstalling the Electrodes and the Insulation Plate

- 1 Reinstall one electrode in the outlet electrode position on the insulation plate.
- 2 Thread the electrode into the outlet position on the insulation plate by hand first.
- 3 Tighten the electrode with the 12 mm hexagonal socket screwdriver.
Do not overtighten. Do not bend the electrodes.
- 4 Holding the insulation plate by the cassette holder, slide the insulation plate under the detector (1.), then tilt it up into place (2.) and push it completely to the left (3.).

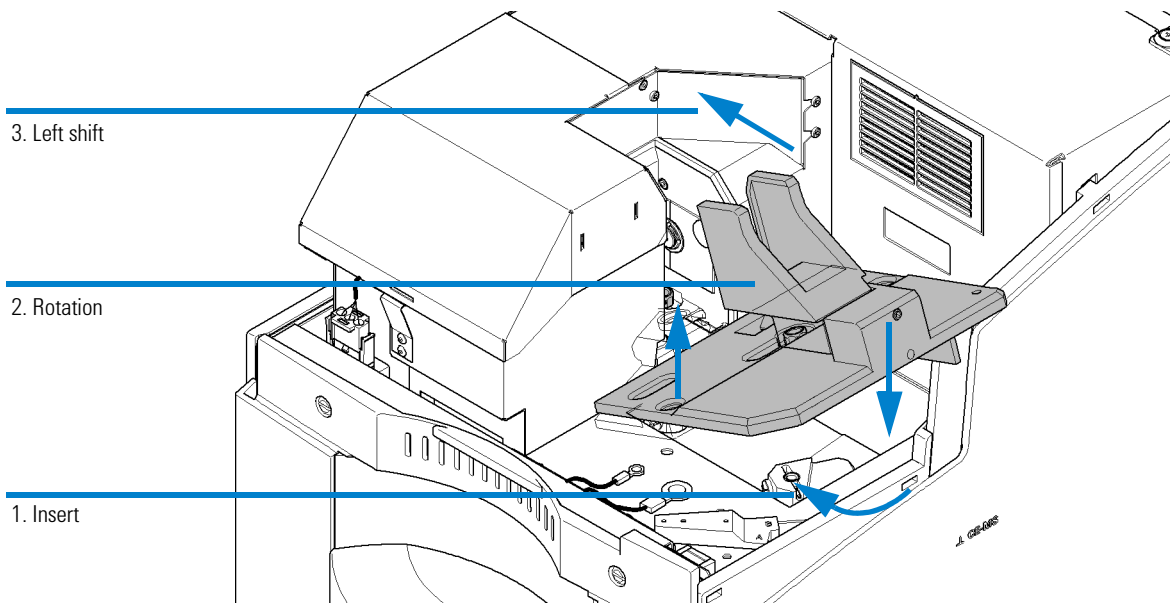
Figure 71 Installation of the insulation plate

Cleaning the Electrodes, Pre-punchers and Insulation Plate

- 5** The insulation plate should be laying flat in position.

The insulation plate is correctly positioned under the detector if the top part of the outlet electrode fits through an oval cut-out.

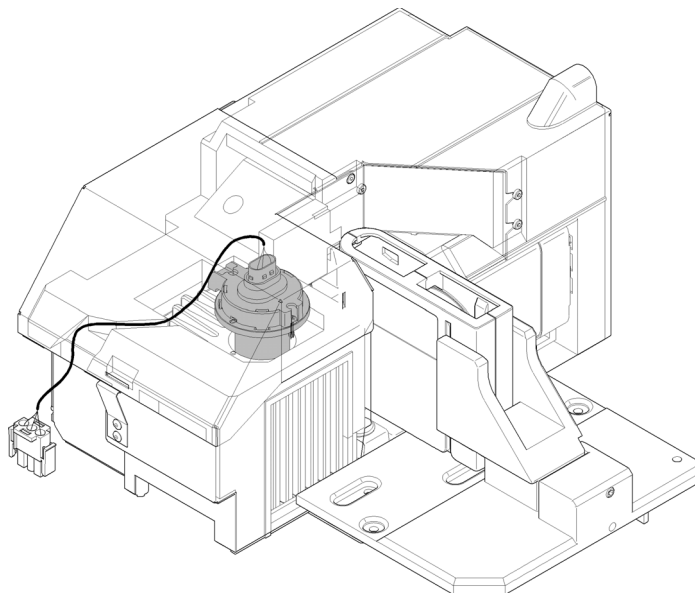
Figure 72 Alignment of the detector

- 6** Secure the insulation plate by reinstalling the two slotted metal screws.
- 7** Reconnect the blue ground wire and fasten the metal screw.
- 8** Insert the inlet electrode into the ring connector of the red high voltage wire. Thread the inlet electrode (and the ring connector) into the insulation plate by hand, then tighten the inlet electrode with the 12 mm hexagonal socket screwdriver.
Do not overtighten.

8 Maintenance

Cleaning the Electrodes, Pre-punchers and Insulation Plate

- 9 Push the liquid handling module back into the instrument as shown in [Figure 74](#) on page 233. To release it from the front position, slightly more force is needed. Carefully move the liquid handling module completely into a position that allows fixation. Please note the metal nose from the liquid handling module which has to fit in a groove of the back metal sheet as shown in [Figure 73](#) on page 232.

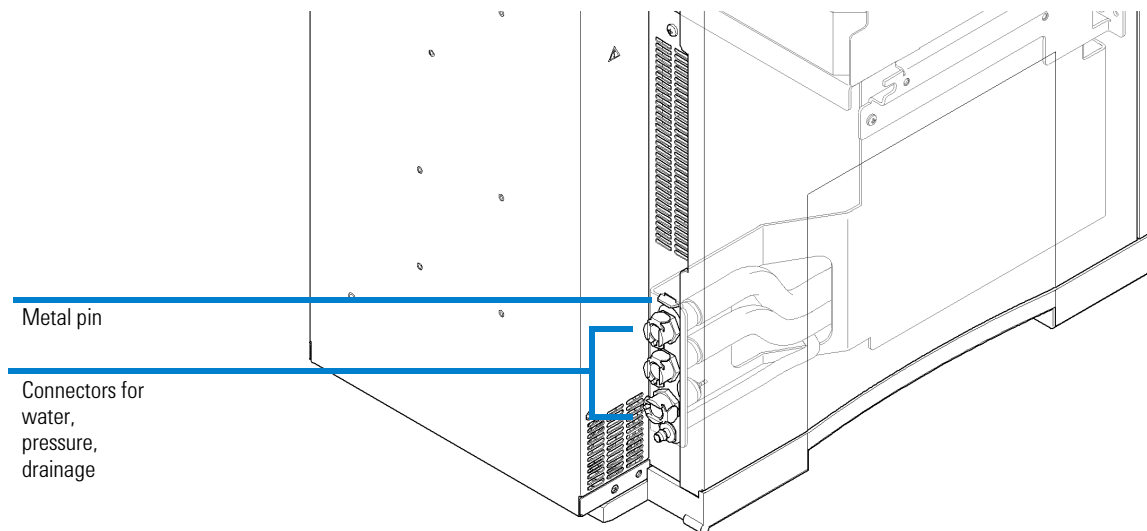
Figure 73 Rear side of liquid handling module

NOTE

Take care not to damage tubes and flat ribbon cables on the right-hand side of the liquid handling module.

Take care that the connectors on the left-hand side, for example for the drainage, fit into their cut-out in the back.

10 Tighten the slotted screws in the front of the liquid handling module.

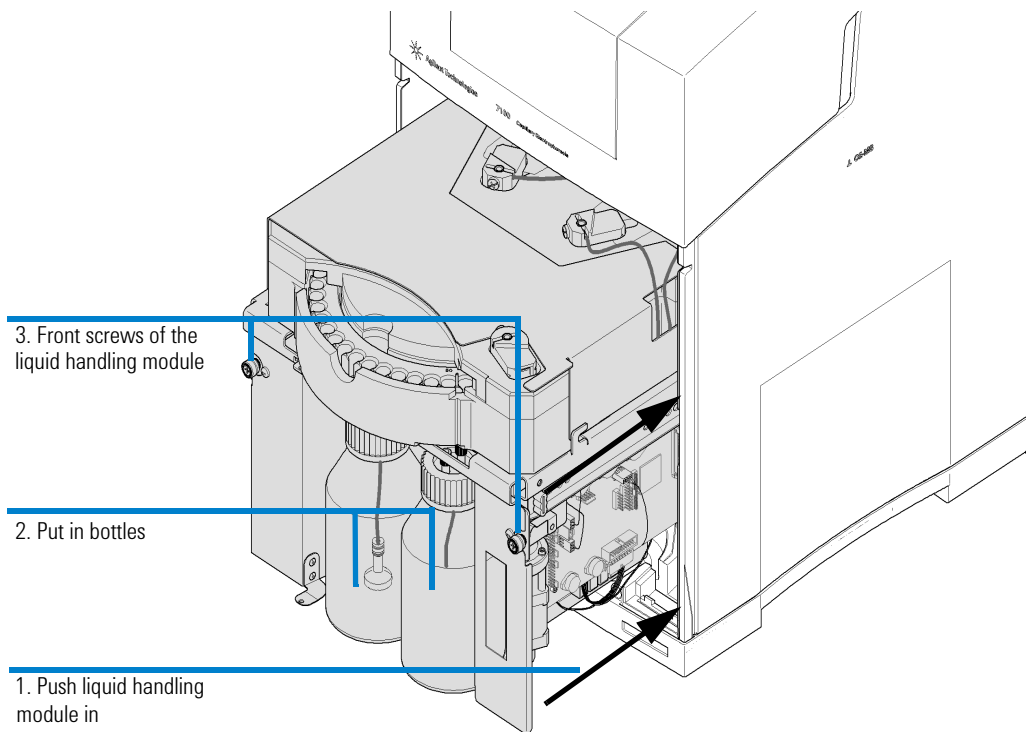
Figure 74 Liquid handling module

11 Reconnect the line power, drainage, external pressure supply, the connectors for an external cooling bath (if applicable), and switch on the instrument.

12 Reestablish the connection to the instrument from the Lab Advisor software. Select **Start monitoring**.

NOTE

No high voltage can be applied, if the front cover is removed. Touching the electrodes inside the instrument is safe. However, under the following circumstances (usage of the alignment center), slow lifter movements with a low force are possible with the following alignment procedure. When you start any lifter movement your hands should be outside the instrument.

13 Open the **CE Control** screen from the **Service & Diagnostic** context > **G7150A CE > CE Control Screen** in the Lab Advisor software. Select the **Alignment Center**. The alignment procedure can be started with the front cover un-installed.

14 Follow the procedure to align the inlet and outlet electrode and the replenishment needle.

The procedure consists of a prepare step and an alignment check.

The alignment is done while the pre-puncher of the respective lift is raised until it is near the electrode or needle tip (prepare-button). Now you can insert the capillary of the electrode alignment tool (**G7100-60107**) from the funnel of the electrode. With careful puncture you can check if the electrode exit points fit exactly to the pre-puncher funnel exit. The electrode or needle might need careful manual bending to make it point towards the exact center of the pre-puncher funnel.

During the alignment check (check-button), the electrode or replenishment needle passes through the funnel several times. An improperly aligned electrode slides along the pre-puncher funnel. The orientation is optimum if the electrode does not move sideways during this dynamic check.

NOTE

Although electrodes will most likely not need alignment thanks to their design, you should still perform the check alignment procedure.

NOTE

The electrodes are fragile. Manual bending during alignment should be done only with great care. Electrodes must be kept straight. Avoid curving.

Reinstall the Capillary Cassette

See [“Inserting the Cassette”](#) on page 145.

Reinstall the Front Cover

- 1 Make sure the front cover screws have a horizontal orientation. They can only click into place in this orientation.
- 2 Keep the front door of the cabinet slightly open.
- 3 Start mounting the front cover by inserting it at the lower end first.
- 4 Carefully push the front cover into the main frame.
You might need to pay special care on the right side due to a connector.
- 5 Close the lid of the instrument.

Maintenance of the Replenishment System

The replenishment system automatically empties and refills the buffer vials in the tray with new buffer from the electrolyte bottle. Typically, replenishment is done before every run or after every 3 to 4 runs, depending on the stability and volume of the running buffer. Initial preparation of the replenishment system is described in [“Preparing the Replenishment System if Needed”](#) on page 125. Please see the general instructions on operations for replenishment maintenance such as flush tubes, clean tubes and clean level sensor ([“Operations for Replenishment Maintenance”](#) on page 239).

Frequency

The replenishment system (bottles and tubing) should be cleaned if:

- using it for the first time,
- changing the buffer, or
- the replenishment system will be idle for some time.

Requirements

- Vials and caps
- Water and/or buffer and isopropanol (depending on the protocol selected)
- ChemStation, Lab Advisor or Instrument Utilities software

Changing the Buffer Composition in the Replenishment System

Please follow these steps.

- 1 Open the Lab Advisor software, connect to the instrument and start monitoring the respective instrument.
- 2 Open the **CE Control** screen from the **Service & Diagnostic** context > **G7150A CE** > **CE Control Screen** and click on **Change Bottles** to release the pressure before you open the bottles.
- 3 Take out the electrolyte bottle and unscrew the cap. The electrolyte bottle has a frit and tubing inside. The frit must not get dirty.
- 4 Fill a minimum 300 mL of buffer into the electrolyte bottle. Make sure the buffer is filtered by using a 0.25 μm or 0.45 μm filter depending on its viscosity.
- 5 Make sure the waste bottle is empty.
- 6 Cap the bottles tightly.
- 7 Choose **Done** in the **Change Bottles** dialog box. Wait until the pressure has built up again.
- 8 To flush the internal tubing system from the electrolyte bottle to the waste bottle including all valves on this pathway, select **Flush Tubes**. During this step the replenishment needle itself is not flushed. Set a flushing time of 0.3 minutes.
- 9 To clean the internal tubing system, including the replenishment, select **Clean Tubes**. During this step the replenishment is used to fill and empty a vial several times. Set the number of repetitions to 5.
- 10 When requested to do so, place a capped and empty vial into a tray position.

The involved volume of the replenishment system is about 5 mL. A 10-time flush is sufficient. The flush time depends primarily on the viscosity of your buffer. Therefore, choose a short time and repeat flush tubes if necessary, otherwise it is possible that the buffer is flushed completely into the waste bottle. This protocol suggests a combination of flush tubes first and a subsequent clean tubes step, with comparably low volume consumption, for a high exchange efficiency and low buffer consumption.

The Replenishment System Will Not Be Used for Some Time

The replenishment system should be flushed first with water and then with isopropanol.

Please follow these steps.

- 1 Open the Lab Advisor software, connect to the instrument and start monitoring the respective instrument.
- 2 Open the **CE Control** screen from the **Service & Diagnostic** context > **G7150A CE > CE Control Screen** and click on **Change Bottles** to release the pressure before you open the bottles..
- 3 Take out the electrolyte bottle and unscrew the cap. The electrolyte bottle has a frit and tubing inside. The frit must not get dirty.
- 4 Fill a minimum 300 mL of water into the electrolyte bottle. Make sure the water is filtered by using a 0.25 µm or 0.45 µm filter depending on its viscosity.
- 5 Make sure the waste bottle is empty.
- 6 Cap the bottles tightly.
- 7 Choose **Done** in the **Change Bottles** dialog box. Wait until the pressure has built up again.
- 8 To flush the internal tubing system from the electrolyte bottle to the waste bottle, including all valves on this pathway, select **Flush Tubes**. During this step the replenishment needle itself is not flushed. Set a flushing time of 2 minutes.
- 9 To clean the internal tubing system, including the replenishment, select **Clean Tubes**. During this step the replenishment is used to fill and empty a vial several times. Set the number of repetitions to 5.
- 10 When requested to do so, place a capped and empty vial into a tray position.
- 11 Execute the clean level sensing operation to clean the replenishment needle entirely.
- 12 Repeat all steps starting with number 2 and use isopropanol instead of water. Empty the waste bottle whenever you fill or refill the electrolyte bottle.

Further Action

If the instrument cannot build-up pressure or a vacuum after performing the above steps, check the following:

- 1 Proper seal of the electrolyte bottle
- 2 Proper seal of the waste bottle
- 3 Tight fittings

If none of these reasons apply, call Agilent Technologies.

Operations for Replenishment Maintenance

The following operations can be started from the ChemStation, Lab Advisor or Instrument Utilities software. From the ChemStation, they can be done by clicking the replenishment vial. It is a vial located between the replenishment bottles in the CE dialog. From the Lab Advisor or Instrument Utilities software, start monitoring, visit the **Tools** section and the **CE Control** screen.

Flush Tubes

The **Flush Tubes** operation directly flushes water, buffer or other cleaning solutions from the electrolyte bottle to the waste bottle. During this task, a reduced pressure is utilized in the electrolyte bottle. The user can set a time for this operation.

The volume in the tubing system from the electrolyte bottle and the waste bottle is about 5 mL. It is recommended that you flush at least with 10-fold volume of fresh buffer. The exchange flow rate of water is approximately 120 to 150 mL/min. Please note the individual viscosity of your buffer system.

It is important to make sure that there is enough buffer (for example a minimum of 300 mL) in the electrolyte bottle for the **Flush Tubes** operation. An excessive **Flush Tubes** time will drain the electrolyte bottle, causing air bubbles to be introduced into the system.

The **Flush Tubes** command does not include a cleaning of the replenishment needle. It requires an execution of the **Clean Tubes** function to clean the last part to the fill needle of the replenishment needle.

Clean Tubes

The **Clean Tubes** operation is meant to replace liquid in tubing from the electrolyte bottle to the replenishment needle and from the replenishment needle to the waste bottle. A vial is lifted to the replenishment needle and filled several times. The number of filling/emptying cycles can be set by users using the Lab Advisor/Instrument Utilities software. This operation is meant for buffer change purposes or for cleaning with water.

It is important to make sure there is enough buffer in the electrolyte bottle for the **Clean Tubes** number of filling/emptying cycles specified (for example a minimum of 50 mL). An excessive **Clean Tubes** time will drain the electrolyte bottle, causing air bubbles to be introduced into the system.

The volume in the tubing system from electrolyte bottle and the waste bottle is approximately 5 mL. Please note that a relatively high amount of buffer is located in the glass frit of the electrolyte bottle. It is recommended that you use the **Flush Tubes** command first to flush the direct tubing connection from the electrolyte bottle to the waste bottle with at least for example 10-times the volume of fresh buffer. Please note the individual viscosity of your buffer system.

NOTE

If the replenishment system will not be used for some time, it should be cleaned first with water and then with isopropanol by the **Flush Tubes** and **Clean Tubes** operations.

Clean Level Sensor

The purpose of the Clean Level Sensor functionality is to dissolve small buffer crystals which may have formed in the tip of the level sensing needle (short needle of the double needle) in the replenishment needle assembly. Remaining pathways of the internal tubing of the level sensing circuit are also cleaned. One indication to use this function is failed level sensing during operation. The other indication is a failed IRP (Injection Replenishment and Pressure) level sensing test from the injection replenishment system tests of the Lab Advisor. You may use this procedure as well if you leave the replenishment system unused for a while.

This function prompts you to load tray position 48 (or user input) with a vial filled with an appropriate cleaning liquid (usually water or current buffer system). When initiated, 8 cleaning cycles are performed.

A cycle consists of:

- 1 Submerging the short needle into the cleaning liquid and briefly sucking cleaning liquid into the short needle.
- 2 Removing the short needle from the cleaning liquid, and sucking a slug of air into the short needle.

The alternate slugs of liquid and air exit the level sensing tubing into the waste bottle. The main screen reply line counts the cycles as they take place. You can check for proper function afterwards with the IRP level sensing test from the injection replenishment system tests or by performing buffer replenishment in a vial.

NOTE

This function is designed to prevent a gradual build-up of buffer crystals. Large, solid plugs of precipitate which have been allowed to form in the short needle may not be removable using this function. In such cases, we recommend exchanging the restrictor capillaries and or the replenishment needle by personnel trained by Agilent Technologies.

NOTE

If the replenishment system will not be used for some time the level sensor should be cleaned with water. Subsequently the replenishment system should be cleaned with water and then with isopropanol by the **Flush Tubes** and **Clean Tubes** operations.

Changing the Air Inlet Filter

Objective

The following procedure is used to replace the air inlet filter. The installed air pump uses ambient air (through this filter) to create pressure inside the electrolyte bottle. The compressed air is needed to flush a capillary and to inject a sample.

Frequency

Air filter replacement should be considered if the environment is very dusty or if the system cannot generate enough pressure. We suggest that you change the air filter at least every three months.

Accessing the Air Filter

The air filter is located inside the electrolyte and waste bottle compartment, see [Figure 75](#) on page 243. Therefore, it is very easy to replace the air filter frequently.

If the instrument cannot build-up either pressure or a vacuum, check the following:

- proper seal of the pressure bottle,
- proper seal of the vacuum bottle, and
- loose fittings on the bottles

If none of these reasons apply, call Agilent Technologies.

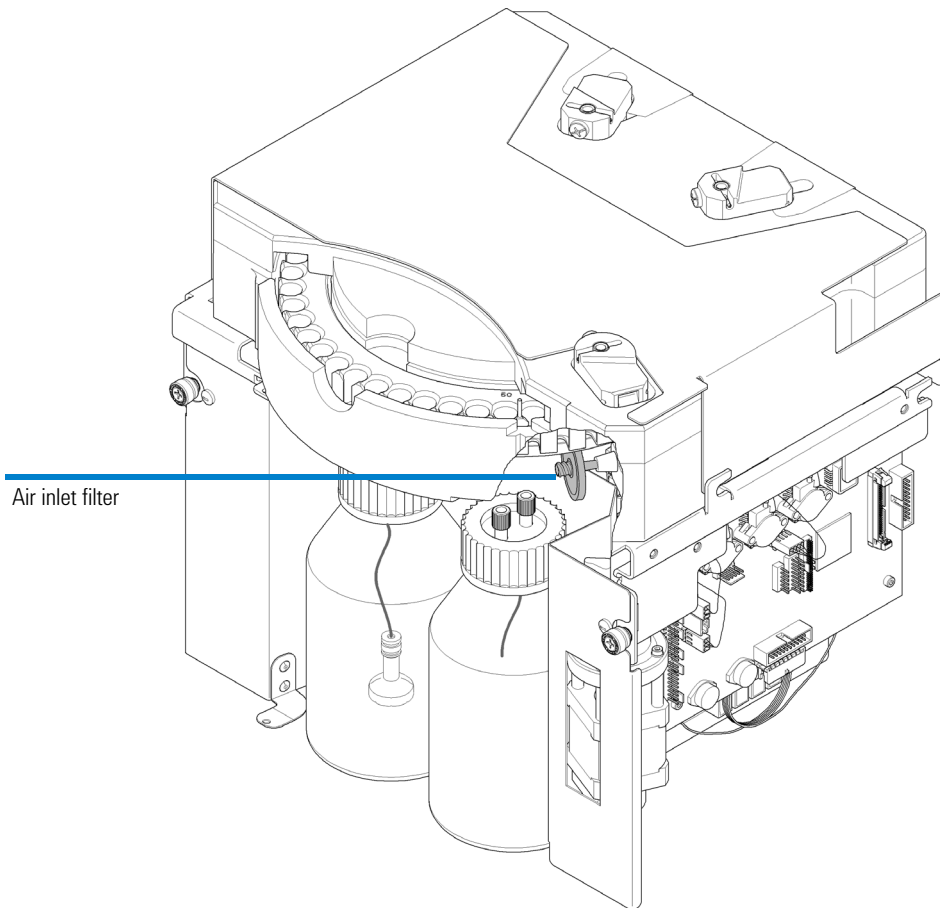
Figure 75 Air inlet filter

NOTE

Do not run the instrument without the air filter installed, since particles may damage the valve system.

Changing the Lamp

The following procedure is used to replace the deuterium lamp in the 7100 CE instrument. Loss of light intensity over time increases baseline noise, which may negatively influence the performance of your analysis.

You should consider exchanging the lamp if a drastic increase in the noise level is encountered which is not caused by any other reason (for example contaminated slit, something absorbing inside the capillary, etc.) or the lamp refuses to ignite. Generally the light intensity drops by 50 % of its original value after approximately 2000 hours of operation. We suggest that you check the lamp intensity before you change the lamp to be sure the noise increase is not due to other reasons (air bubbles inside the capillary, alignment interface or detection window contaminated).

Please use the **Intensity Test** in the **Detector Test** suite in the Lab Advisor/Instrument Utilities software to determine the lamp intensity. Check the Lab Advisor/Instrument Utilities software to read out the lamp tag information on the current lamp lifetime.

It may happen that a hot lamp refuses to ignite the first time. Please wait until the lamp has cooled down.

NOTE

Only lamps provided by Agilent Technologies are supported in the 7100 CE instrument. The lamp tag stores various data and details on the ignition program of the deuterium lamp. A lamp without a tag will fail to ignite and could lead to instrument errors.

Removing the Existing Lamp

Parts required Pozidriv screwdriver (part of the accessory kit).
Deuterium lamp for the Agilent CE instrument.

- 1 Select **Exit** from the **File** menu to exit the Agilent ChemStation. Exit both online and offline copies.
- 2 Turn off the line power to the Agilent CE instrument.
- 3 Open the top cover.
- 4 Open the detector cover.

8 Maintenance

Changing the Lamp

- 5 Disconnect the lamp cable's 8-fold connector. Pull the connector, not the cable.

WARNING

Injuries due to burns

Touching a hot lamp can cause serious burns.

- If you were using the instrument shortly before, the lamp may be very hot. Wait until it has cooled down before touching the lamp.

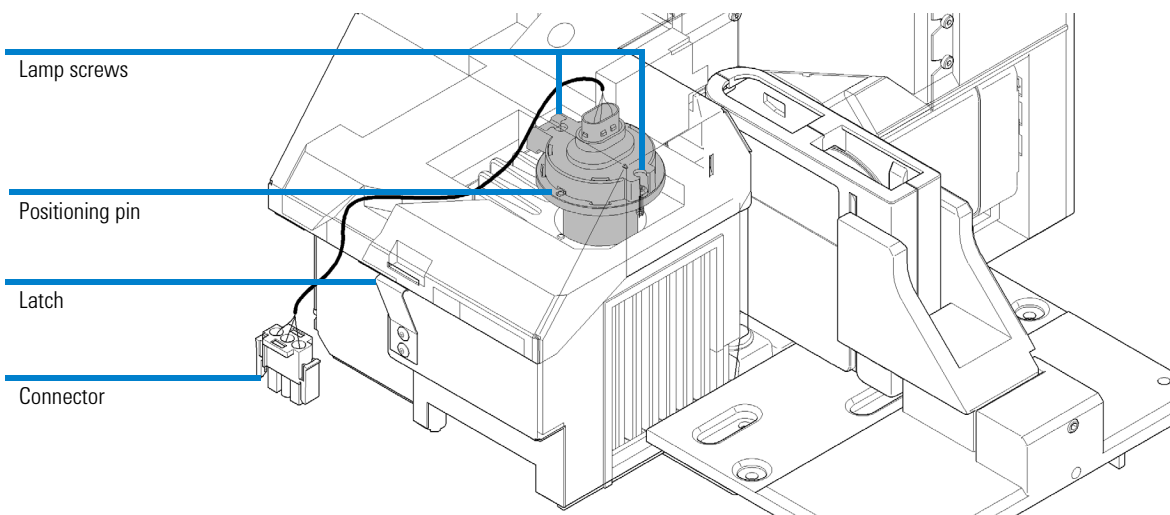
Figure 76 Exchanging a lamp

- 6 Loosen the two lamp screws using the Pozidriv screwdriver and lift the lamp out carefully.

Inserting a New Lamp

- 1 Hold the new lamp by the metal ring to avoid touching the glass surface. Fingerprints absorb light and can reduce light intensity.
- 2 Insert the lamp into the lamp compartment. Position the lamp so that the notch in the lamp ring matches with the pin on the detector housing.
- 3 Tighten the two lamp screws.
- 4 Connect the lamp cable connector and fix the cables into the channel in the foam.
- 5 Close the detector lid without clamping the lamp cable.
- 6 Close the top cover. Turn line power on.
- 7 Boot the Agilent ChemStation and turn the lamp on.
- 8 Allow the lamp to warm up for about 1 hour.
- 9 Perform an **Intensity Test** in the **Detector Test** suite in the Lab Advisor/Instrument Utilities software to determine the initial lamp intensity of your new lamp. Store the result for your own reference.

Further Action

If the instrument fails this test, check the following:

- correct seating and connection of the lamp cables,
- correct seating of the cassette, and
- if the green alignment interface without capillary is used, check for contamination of the interface.

If none of these reasons apply, call Agilent Technologies.

Cleaning the Instrument

The instrument case should be kept clean. Cleaning should be done with a soft cloth slightly dampened with water or a solution of water and mild detergent. Do not use an excessively damp cloth as liquid may drip into the instrument.


WARNING

Risk of instrument damage

Liquid dripping into the electronic compartment of your module.

Liquid in the module electronics can cause shock hazard and damage the module.

- Do not use an excessively damp cloth during cleaning.
 - Drain all solvent lines before opening any fittings.
-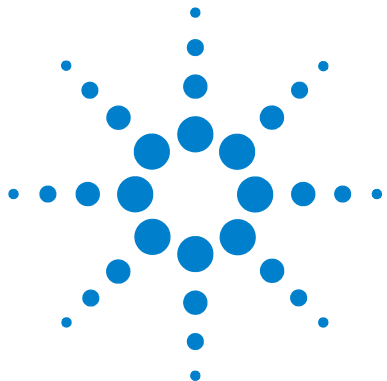

9 Parts and Materials for Maintenance


Overview of Maintenance Parts [250](#)

This chapter provides information on the parts required for maintenance.

Overview of Maintenance Parts

Table 17 Maintenance parts

Description	Part Number
Long life Deuterium lamp (8-pin)	5190-0917
Electrode assembly, standard	G7100-60007
Electrode assembly, short	G7100-60033
Electrode O-ring, silicone 5/pk	5062-8544
Electrolyte bottle, 500 mL, Pressure Plus	9301-6531
Electrolyte bottle, 100 mL, Pressure Plus	9301-6532
Electrolyte bottle cap	9300-1747
Bottle sealing O-ring	0905-1163
Glass filter, solvent inlet	5041-2168
Filter frit adapter, 3 mm, 4/pk	5062-8517
Pre-puncher	G1600-67201
Screws for pre-puncher, 10/pk	G1600-62402
Air filter, 5 µm	3150-0619
Capillary cassette	G7100-60002
Optical filter for DAD, 260 nm	G7100-62700
Optical filter for DAD, 280 nm	G7100-68750
Vials	See Table 13
Alignment interfaces	See Table 14
Fuses	See “ 7100 Capillary Electrophoresis Instrument Electronics ” on page 193

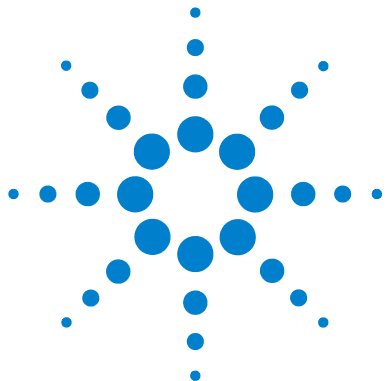10 Appendix

General Safety Information	252
Allgemeine Sicherheitsinformation	255
Informations générales de sécurité	258
Informazioni generali sulla sicurezza	261
Información de seguridad	264
安全に関する一般的な情報	267
一般安全信息	270
Setting Up a Test Sample Run	272
Chemical and Biological Safety	276
The Waste Electrical and Electronic Equipment (WEEE) Directive (2002/96/EC)	277
Radio Interference	278
Sound Emission	279
Legal Notice	280
Agilent Technologies on Internet	281

This chapter provides additional information regarding safety, legal considerations and Agilent information on the Internet.

This is a Safety Class I instrument (provided with terminal for protective earthing) and has been manufactured and tested according to international safety standards.

General Safety Information

The following general safety precautions must be observed during all phases of operation, service, and repair of this instrument. Failure to comply with these precautions or with specific warnings elsewhere in this manual violates safety standards of design, manufacture, and intended use of the instrument. Agilent Technologies assumes no liability for the customer's failure to comply with these requirements.

General

This is a Safety Class I instrument (provided with terminal for protective earthing) and has been manufactured and tested according to international safety standards.

Operation

Before applying power, comply with the installation section. Additionally the following must be observed.

Do not remove instrument covers when operating. Before the instrument is switched on, all protective earth terminals, extension cords, auto-transformers, and devices connected to it must be connected to a protective earth via a ground socket. Any interruption of the protective earth grounding will cause a potential shock hazard that could result in serious personal injury. Whenever it is likely that the protection has been impaired, the instrument must be made inoperative and be secured against any intended operation.

Make sure that only fuses with the required rated current and of the specified type (normal blow, time delay, and so on) are used for replacement. The use of repaired fuses and the short-circuiting of fuse holders must be avoided.

Some adjustments described in the manual, are made with power supplied to the instrument, and protective covers removed. Energy available at many points may, if contacted, result in personal injury.

Any adjustment, maintenance, and repair of the opened instrument under voltage should be avoided whenever possible. When inevitable, this has to be carried out by a skilled person who is aware of the hazard involved. Do not attempt internal service or adjustment unless another person, capable of rendering first aid and resuscitation, is present. Do not replace components with power cable connected.

Do not operate the instrument in the presence of flammable gases or fumes. Operation of any electrical instrument in such an environment constitutes a definite safety hazard.

Do not install substitute parts or make any unauthorized modification to the instrument.

Capacitors inside the instrument may still be charged, even though the instrument has been disconnected from its source of supply. Dangerous voltages, capable of causing serious personal injury, are present in this instrument. Use extreme caution when handling, testing and adjusting.

When working with solvents please observe appropriate safety procedures (e.g. safety glasses, gloves and protective clothing) as described in the material handling and safety data sheet by the solvent vendor, especially when toxic or hazardous solvents are used.

Safety Symbols

Table 18 Safety Symbols

Symbol	Description
	The apparatus is marked with this symbol when the user should refer to the instruction manual in order to protect risk of harm to the operator and to protect the apparatus against damage.
	Indicates dangerous voltages.
	Indicates a protected ground terminal.
	Indicates eye damage may result from directly viewing the light produced by the deuterium lamp used in this product.
	The apparatus is marked with this symbol when hot surfaces are available and the user should not touch it when heated up.

WARNING

A WARNING

alerts you to situations that could cause physical injury or death.

- Do not proceed beyond a warning until you have fully understood and met the indicated conditions.

CAUTION

A CAUTION

alerts you to situations that could cause loss of data, or damage of equipment.

- Do not proceed beyond a caution until you have fully understood and met the indicated conditions

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Setting Up a Test Sample Run

This instrument check demonstrates that the instrument is correctly installed and connected. Please use a fused silica capillary with 40 cm effective length and 50 µm inner diameter from the HPCE start-up capillary kit (part number **G1600-68700**) with the green alignment interface (part number **G7100-60210**). The instrument should be in CE mode.

Load chemicals from the Installation Qualification Chemical kit (part number **5063-6514**) according to the table below. The Installation Qualification Chemical kit contains:

- Buffer (20 mM borate, pH 9.3, 100 mL)
- Test sample (4-hydroxyacetophenone 1 mM, 3 mL)
- Conditioning solution (1.0 N sodium hydroxide, 100 mL)

Table 25 Solution for Checkout Run

Location	Solution	Volume
Vial 1	1.0 N sodium hydroxide	500 µL
Vial 2	Water	200 µL
Vial 4	Buffer	500 µL
Vial 5	Buffer	500 µL
Vial 6	Buffer	500 µL
Vial 7	Test sample	500 µL
Vial 8	Water/waste	200 µL

Choose the menu item **Method > New method** to prepare a method with the parameters given below. Save the edited method under a new name. Choose the menu item **RunControl > Sample Info...** and enter the operator name, directory and file name. Enter Vial location 7, enter sample name “1 mM pHAP” and click **Run method** to start the analysis.

A typical electropherogram will show a significant peak at a migration time of around 3 minutes.

Method Parameter CE (Default Settings)

Vials

- Inlet home vial 4
- Outlet home vial 5

Cassette Temperature

- 20.0°C

High Voltage System

- Enable high voltage
- Voltage 0.0 V (will ramp up in accordance with timetable)
- Current 300 μ A
- Power 6.0 W
- Low current limit OFF

Stoptime

- 10.0 minutes

Posttime

- OFF

Replenishment

- Replenishment and preconditioning parallel processing
- No replenishment used

Preconditioning

CE Conditioning

- Flush inlet 1, outlet 8 for 5 minutes
- Wait 5 minutes
- Flush inlet 2, outlet 8 for 5 minutes
- Flush inlet 6, outlet 8 for 10 minutes

CE Injection

- Inject by pressure, 50 mbar, 4 seconds

Timetable

- Change voltage 0.2 minutes, 30 kV

Postconditioning

- No postconditioning used

Method Parameter DAD

Signals: Signal, Bw (nm) Reference, Bw

- Use signal A: Yes wavelength 325, Bandwidth 10, no Reference
- Use signal B: Yes wavelength 192, Bandwidth 4, no Reference
- Use signal C: Yes wavelength 235, Bandwidth 10, no Reference

Peak width

- > 0.1 minutes (2.0 s response time) (2.5 Hz)

Stoptime

- As with CE

Posttime

- Off

Spectrum

- None

Autobalance

- Prerun Autobalance

Lamps Required for Acquisition

- UV lamp

Timetable

- No entry

Chemical and Biological Safety

WARNING

Toxic and hazardous solvents

The handling of solvents and reagents can hold health risks.

- When working with solvents observe appropriate safety procedures (for example, goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the solvent vendor, especially when using toxic or hazardous solvents.
-

WARNING

Pathogenic, toxic, radioactive samples and genetically modified organisms

If pathogenic, toxic, or radioactive samples are intended to be used in this instrument, it is the responsibility of the user to ensure that all necessary safety regulations, guidelines, precautions and practices are adhered to accordingly. This includes also the handling of genetically modified organisms. Ask your laboratory safety officer to advise you about the level of containment required for your application and about proper decontamination or sterilization procedures to follow if fluids escape from containers.

- Observe all cautionary information printed on the original solution containers prior to their use.
 - Because leaks, spills, or loss of sample may generate aerosols, observe proper safety precautions.
 - Agilent 7100 CE instrument covers have not been designed as bioseals for aerosols or liquid containment.
 - Handle body fluids with care because they can transmit disease. No known test offers complete assurance that they are free of microorganisms. Some of the most virulent – Hepatitis (B and C) and HIV (I-V) viruses, atypical mycobacteria, and certain systemic fungi – further emphasize the need for aerosol protection.
 - Always follow local state and federal biohazard handling regulations when disposing of biohazardous waste material.
 - Handle all infectious samples according to good laboratory procedures and methods to prevent spread of disease.
 - Dispose of all waste solutions and products according to appropriate environmental, health and safety guidelines.
-

The Waste Electrical and Electronic Equipment (WEEE) Directive (2002/96/EC)

Abstract

The Waste Electrical and Electronic Equipment (WEEE) Directive (2002/96/EC), adopted by EU Commission on 13 February 2003, is introducing producer responsibility on all Electric and Electronic appliances from 13 August 2005.

NOTE

This product complies with the WEEE Directive (2002/96/EC) marking requirements. The affixed label indicates that you must not discard this electrical/electronic product in domestic household waste.

Product Category: With reference to the equipment types in the WEEE Directive Annex I, this product is classed as a "Monitoring and Control instrumentation" product.

Do not dispose off in domestic household waste.

To return unwanted products, contact your local Agilent office, or see www.agilent.com for more information.

Radio Interference

Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.

Test and Measurement

If test and measurement equipment is operated with equipment unshielded cables and/or used for measurements on open set-ups, the user has to assure that under operating conditions the radio interference limits are still met within the premises.

Sound Emission

Manufacturer's Declaration

This statement is provided to comply with the requirements of the German Sound Emission Directive of 18 January 1991.

This product has a sound pressure emission (at the operator position) < 70 dB.

- Sound Pressure $L_p < 70$ dB (A)
- At Operator Position
- Normal Operation
- According to ISO 7779:1988/EN 27779/1991 (Type Test)

Legal Notice

The Agilent 7100 Capillary Electrophoresis System is sold for research only. By purchasing this instrument, the purchaser is granted the limited right to use only this instrument. Purchase of this instrument does not include any right, express or implied, to use any other patented product, method or process, or to use any other portion or component of any patented system or systems, software, microfluidic devices or reagents, either alone or conjunction with this product, unless use of such method, process or other portion or component is separately authorized.

Agilent Technologies on Internet

For the latest information on products and services visit our worldwide web site at:

<http://www.agilent.com>

Select **Products/Chemical Analysis**

It will provide also the latest firmware of the Agilent 7100 Capillary Electrophoresis System for download.

10 Appendix
Agilent Technologies on Internet

Index

A

accessing, 219
accessories, 69, 250
acquisition
 current, 119
 detector signals, 121
 power, 113
 pressure, 119
 spectra, 122
 temperature, 111
 voltage, 112
AD converter, 201, 206
alignment interface, 137, 211
 color coding, 137
 installation of capillary, 137
 removing, 135, 145
autobalance, 123

B

buffer, 125
 buffer vials, 129
 change, 125
 change during run, 125
 composition, 237
 liquid level, 131
 preparation, 125
 replenishment, 125

C

capillary
 installation into cassette, 141
capillary cassette
 opening, 141
 removing, 135
 temperature, 111
capillary conditioning, 115

capillary detection window, 138
capillary electrochromatography, 169
capillary electrophoresis test kit, 272
capillary insertion tool, 138
capillary lock, 142
carousel, 143
cassette, 142
 closing, 144
 installation, 145
 installation of capillary, 142
 removing from instrument, 135
 temperature, 111
cassette lock, 142
CE diagram, 98
CE timetable, 119
change cassette function, 135
changing, 246
changing of fuses, 202
checkout run, 272
ChemStation, 93
 starting, 95
cleaning, 211, 216, 225
condensing water, 20, 111
conditioning, 115
corresponding alignment interface
 detection window, 138
 extended light path, 137
 installation into alignment
 interface, 138
corresponding alignment interfaces, 137
corrosion, 126
current, 112, 119

D

detector, 121
 diagram, 100
 icon, 100
 lamp, 100, 250
 parameters, 108, 121
 posttime, 122
 signals, 123
 stoptime, 122
 timetable, 123
detector cover, 192
diagnostic, 189

E

electric parameters, 112
electrodes, 112, 225
electrolyte bottle, 101
environment, 18, 20
 condition, 20
extended light path capillaries, 138

F

firmware, 196, 197
flush tubes, 239
fraction collection, 148
front cover, 218

G

gradients, 119

H

High Pressure
 home vials, 129
 humidity, 111

Index

high Pressure
 home vials, 110
high pressure, 157, 169
humidity, 20

I

inlet home vial, 129
installation, 230
instrument, 92, 101
 starting, 93
instrument control, 93
insulation plate, 227

L

Lab Advisor, 189
lamp
 switching on, 124
leaks, 183
levelsensor, 240
line frequency, 21
line fuse holder, 193
line power, 72
line voltage, 21
liquid level, 129
 buffer vials, 132
 outlet vials, 133
 sample vials, 131
 waste vials, 133
lower alarm limit, 113

M

main system, 196
maintenance, 209
 alignment interface, 137
method parameters, 106, 273

N

noise level, 121, 279

P

paragraph tags, 18, 68, 92, 192, 210, 250, 251
peak width, 121
polarity, switching, 112
posttime, 114
power, 113
power supply, 21
pre-puncher, 226
 cleaning, 226
 installation, 223
prepuncher
 installation, 223
pressure, 117

R

raw data, 196
reference wavelength, 121, 275
removing, 245
removing cassette, 135
replenishment system, 114
 installation of bottles, 125
 preparation, 125
resident system, 196

S

sample vials, 131
 liquid level, 131
 preparation, 131
SDS, 188
setup CE method, 106
signals, 121
spectra, 122
starting instrument, 95, 105
stoptime, 114
straight capillaries, 137
switching during run, 123

T

tags
 paragraphs, 18, 68, 92, 192, 210, 250, 251
Telnet, 86
temperature, 111
 environment, 111
test
 instrument, 189
timetable, 119
tray, 101
tray cooling, 204
 manual operation, 134
tubes
 clean, 240

U

unpacking, 68
updates, 197

V

vial sensor, 134
vials, 129
voltage, 21, 112

W

waste bottle, 128
waste vials, 133
water bath, 204
wavelength, 121
wavelength calibration, 147

www.agilent.com

In This Book

This manual contains technical reference information about the Agilent 7100 Capillary Electrophoresis System.

The manual describes the following:

- introduction and specifications,
- installation,
- using and optimizing,
- troubleshooting,
- maintenance
- parts identification,
- safety and related information.

© Agilent Technologies 2009 - 2019

Printed in Germany
11/2019

G7100-90000

Rev. C.

SD-29000166 Rev. C

Agilent Technologies