VWR® Ministar & MiniStar blueline Centrifuge

INSTRUCTION MANUAL

European Catalogue Numbers:

VWR MiniStar

Euro Plug: 521-2161

UK Plug: 521-2162

Swiss Plug: 521-2163

VWR MiniStar blueline

Euro Plug: 521-2319

UK Plug: 521-2320

Swiss Plug: 521-2321

Version: 2

Issued: 17, 08, 2017

Legal Address of Manufacturer

Europe

VWR International bvba

Researchpark Haasrode 2020

Geldenaaksebaan 464

B-3001 Leuven

+ 32 16 385011

<http://be.vwr.com>

Country of origin: Korea

Table of Contents

Symbols and Conventions.....	1
Safety Information	1
Package Contents	1
Installation	2
Intended Use	2
Product Specifications.....	2
Overview.....	2
Operation.....	3
Repair and Maintenance	4
Accessories and Spare Parts	4
Technical service	5
Warranty	5
CE Certificate	6
Equipment Disposal.....	8

Symbols and Conventions

The following chart is an illustrated glossary of the symbols that are used in this manual

	<p>CAUTION This symbol indicates a potential risk and alerts you to proceed with caution.</p>
	<p>CAUTION This symbol indicates the presence of high voltage and warns the user to proceed with caution.</p>
	<p>This symbol identifies a Protective Earth (PE) terminal, which is provided for connection of the supply system's protective earth (green or green/yellow) conductor.</p>

Safety Information

NEVER use the centrifuge in any manner not specified in these instructions.

NEVER operate the centrifuge without a rotor properly attached to the shaft

NEVER fill tubes while they are in the rotor. Liquid spillage may harm unit.

NEVER put hands in the rotor area unless the rotor is completely stopped.

NEVER move the centrifuge while the rotor is spinning.

NEVER use solvents or flammables near this or other electrical equipment.

NEVER centrifuge flammable, explosive or corrosive materials

NEVER centrifuge hazardous materials outside of a hood or proper containment facility

ALWAYS load the rotor symmetrically. Each tube should be counterbalanced by another tube.

Weight of individual tubes may not exceed 3.0g.

ALWAYS locate the centrifuge within easy access to an electrical outlet.

ALWAYS use only microcentrifuge tubes made from plastic and designed to withstand centrifugal forces of at least 2,000 xg

Please read this manual in its entirety to ensure safe operation of the unit.

Package Contents

Description	Quantity
8 Place tube rotor	1 each
4 place PCR strip rotor	1 each
0.5mL tube adapters	Pack of 8
0.4mL tube adapters	Pack of 8
Power cord	1 each
User Manual	1 each

Installation

Bench top area required for the clearance envelope for the intended use

Intended Use

- For Quick Spin-down from Walls or Caps of Centrifuge Tubes
- Complete Set with Circular Fixed-Angle Rotor and Strip Rotor

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

Product Specifications

Dimensions	13.2 x 13.2 x 11.2 cm
Weight	0.55 kg
Electrical	AC 230V, 50.60Hz, 65mA
Fuse	0.5A, 250V (F)
Speed	6000 rpm
RCF	2000 x g
Acceleration time in s	5
Deceleration time in s	1.5
Noise level in db (A)	56
Power consumption in W	5

Overview

The VWR MiniStar Centrifuge is supplied with two rotors, one for microtubes and one for 0.2ml PCR strip tubes. The 8 place rotor is designed to centrifuge up to eight individual 1.5/2.0ml plastic microcentrifuge tubes. It will also accept 0.5ml and 0.4ml tubes with the adapters supplied with the unit. Adapters are available separately for 0.2ml tubes. The strip tube rotor is designed to centrifuge four 8 x 0.2ml strip tubes. It will also accept up to 32 individual 0.2ml tubes. Both rotors are designed for applications requiring relatively low g-forces, such as microfiltration, cell separation and quick spin downs from the walls of tubes.

Please read this manual thoroughly before attempting to operate the centrifuge.

Operation

Before operating the VWR MiniStar centrifuge, be sure that the power switch on the back of the unit is in the “on” position. To begin a run, simply close the lid of the centrifuge. No other controls are required. To stop rotation, press and release the tab on the front of the lid. This will release the lid lock, allowing the lid to self open and the rotor to come to a gradual stop. Caution: Do not attempt to open the lid or remove samples until the unit has come to a complete stop. Once the lid has fully opened and the rotor has stopped, samples may be removed from the rotor.

Safe operation of the centrifuge requires that the rotor be loaded in a balanced fashion. Failure to load the rotor properly presents a hazardous condition and may damage the centrifuge. Individual tubes, strip tubes and/or adapters must always be loaded symmetrically to ensure proper balance. Never run the centrifuge with only one strip or tube in place.

The VWR MiniStar centrifuge is supplied with two interchangeable quick release rotors. No tools are required to exchange the rotors or attach them to the motor shaft. The rotors can be easily exchanged with one hand by way of the quick release system. Removal of the rotors is done by placing your thumb on top of the motor shaft and lifting up with two fingers under the rotor. Place the other rotor into position on the motor shaft and snap in.

(8-place tube rotor and 4-place strip rotor)

Repair and Maintenance

Instructions specify inspection of any means of fixing the equipment to the mounting surface and the condition of the mounting surface itself

Instructions specify safeguards for the operator during cleaning

Inspection of the protective casing

Frequency of inspection and the method of replacement of BIOSEALS and other bio-containment components

The instructions for use state the precautions when the materials to be used with a laboratory centrifuge are known to be toxic, radioactive, or contaminated with pathogenic micro-organisms

Before using any cleaning or decontamination methods except those recommended by the manufacturer, check with the manufacturer that the proposed method will not damage the equipment. To clean the centrifuge, use a damp cloth and a mild, noncorrosive detergent (pH <8).

After cleaning, ensure all parts are dried thoroughly before attempting to operate the unit. Do not immerse the centrifuge in liquid or pour liquids over it.

Accessories and Spare Parts

Description	Quantity	Cat. No.
PCR Strip Rotor, 4-places	1	521-2210 (blue)
		521-2322 (white)
Tube Rotor, 8-places	1	521-2211 (blue)
		521-2323 (white)
Adapter 0,5mL tubes	Pack of 6	521-2817
Adapter 0,4mL	Pack of 6	521-2818

Technical service

Web Resources

Visit the **VWR** website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit.

www.vwr.com.

Warranty

VWR warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, **VWR** will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the product in its local environment. **VWR** will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

VWR International bvba
Haasrode Researchpark 2020, Geldenaaksebaan 464, B-3001 Leuven
<http://be.vwr.com>

EU Declaration of Conformity

Manufacturer: VWR International bvba, Haasrode Researchpark 2020, Geldenaaksebaan 464, B-3001 Leuven

The undersigned states that this declaration is issued under the sole responsibility of the manufacturer, VWR International bvba, and that the object(s) of the declaration described in annex 1 are in conformity with the relevant Union harmonization legislation.

In the event of unauthorized modification of any of the products listed in the annex 1, this declaration becomes invalid.

Object Name: VWR MiniStar Centifuge

Relevant EC Directives:

Low Voltage Directive	2014/35/EU
EMC Directive	2014/30/EU
RoHS	2011/65/EU

Harmonized Standards:

EN 61326-2-1:2013
EN 61010-1:2010
EN 61010-2-020:2013

Document Number: 21264.2017.01

Place & date of issue: Leuven, 28/08/2017

Signed for and on behalf of VWR International bvba:

Yves van Damme
European Regulatory Affairs Manager

Kris Nijs
Managing Director

Tel: 016 385 011 – Fax: 016 385 385 Email : vwr.be@vwr.com
BTW.TVA: BE0403593.343

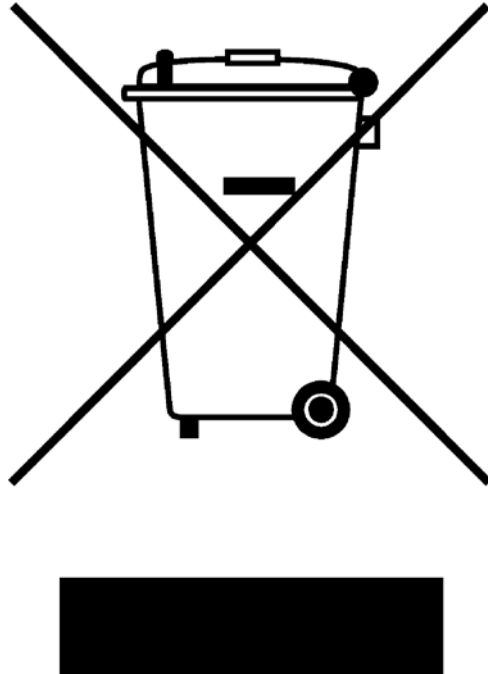
Annex 1

Cat. No.	Description
521-2161	VWR MiniStar, EU-Plug
521-2162	VWR MiniStar, UK-Plug
521-2163	VWR MiniStar, CH-Plug
521-2319	VWR MiniStar BlueLine, EU-Plug
521-2320	VWR MiniStar BlueLine, UK-Plug
521-2321	VWR MiniStar BlueLine, CH-Plug

M3-F46_V2

Equipment Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
Email: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR
International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email sales.ie@vwr.com

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Teknolojileri Ltd. Şti.
VWR International China Co., Ltd
VWR Singapore Pte Ltd